
JE

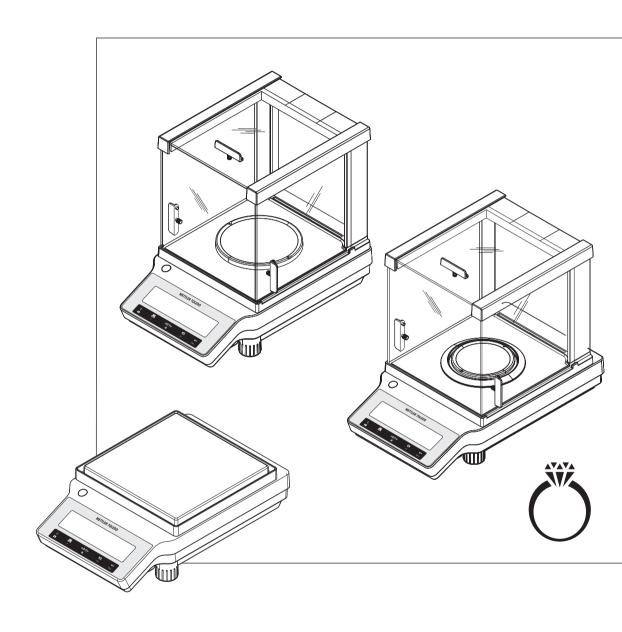



# **Table of Contents**

1	Introdu 1.1	ction Conventions and symbols used in these Operating Instructions	<b>3</b>
2	<b>Safety</b> 2.1 2.2	Information Definition of signal warnings and symbols	<b>4</b> 4
3	<b>Desigr</b> 3.1	Overview	7 7 8 9
4	Install 4.1 4.2 4.3 4.4 4.5 4.6	Unpacking and delivery inspection Installing components Installing protective cover Selecting a location Connecting the balance Setting up the balance 4.6.1 Switching on the balance 4.6.2 Leveling the balance 4.6.3 Setting date and time 4.6.4 Adjusting the balance Adjustment (calibration) 4.7.1 Adjustment with internal weight 4.7.2 Adjustment with external weight 4.7.3 Customer fine adjustment	12 13 15 16 16 17 17 17 17 19 19 20 21
	4.8 4.9 4.10	Performing a simple weighing Weighing below the balance Transporting the balance	22 23 24
5	The Me 5.1 5.2	What is in the menu?  Description of menu topics.  5.2.1 Main menu  5.2.2 Basic menu  5.2.3 Advanced menu  5.2.4 Interface menu	25 26 26 26 27 29
6	<b>Applic</b> 6.1 6.2 6.3 6.4 6.5 6.6 6.7 6.8	Application piece counting Application percent weighing Application check weighing Application statistics Application totaling Application multiplication factor weighing Application division factor weighing Application division factor weighing Application density 6.8.1 Density determination of solids 6.8.2 Density determination of liquids 6.8.3 Formulae used to calculate density	344 377 399 411 433 444 466 488 500 500
7		unication with Peripheral Devices	<b>53</b>

8	Error	and Status Messages	55
	8.1	Error messages	55
	8.2	Status messages	56
9	Maint	enance	57
	9.1	Cleaning and service	57
	9.2	Cleaning the draft shield	58
	9.3	Disposal	58
	9.4	Firmware (Software) Updates	58
10	Techn	ical Data	59
	10.1	General data	59
	10.2		61
		10.2.1 Carat balances with readability of 0.001 ct / 0.1 mg	
		10.2.2 Gold balances with readability of 1 mg	
		10.2.3 Gold balances with readability of 10 mg	
	10.3	Dimensions	
		10.3.1 Carat balances with readability of 0.001 ct / 0.1 mg	
		10.3.2 Gold balances with readability of 1 mg	
		10.3.3 Gold balances with readability of 10 mg	
	10.4		
		10.4.1 RS232C interface	
		10.4.2 MT-SICS interface commands and functions	
11	Acces	sories and Spare Parts	72
	11.1	Accessories	72
	11.2		76
	Index		79

Table of Contents Jewelry Balances

# 1 Introduction

Thank you for choosing a METTLER TOLEDO balance. The balances of the Jewelry line combine a large number of weighing possibilities with easy operation.

The operating instructions are based on the initially installed terminal firmware (software) version V 1.01.

www.mt.com/jewelry

3

# 1.1 Conventions and symbols used in these Operating Instructions

Key and / or button designations are shown in graphic or bold text (e.g.  $\blacksquare$ ). These symbols indicate an instruction:

- prerequisites
- 1 steps
- 2 ...
- ⇒ results



This symbol indicates press key briefly (less than 1.5 s).



This symbol indicates press and hold key down (longer than 1.5 s).



This symbol indicates a flashing display.

Jewelry Balances Introduction

# 2 Safety Information

- Read and understand the instructions in this manual before you use the balance.
- Keep this manual for future reference.
- Include this manual if you pass on the balance to other parties.

If the balance is not used according to the instructions in this manual or if it is modified, the safety of the user may be impaired and Mettler-Toledo GmbH assumes no liability.

### 2.1 Definition of signal warnings and symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

**WARNING** for a hazardous situation with medium risk, possibly resulting in death or severe

injury if not avoided.

**CAUTION** for a hazardous situation with low risk, resulting in minor or moderate injury if not

avoided.

**NOTICE** for a hazardous situation with low risk, resulting in damage to the balance, other

material damage, malfunctions and erroneous results, or loss of data.

Note (no symbol)

for useful information about the product.



General hazard



Electrical shock



NOTICE

# 2.2 Product safety information

Your balance represents state-of-the-art technology and complies with all recognized safety rules, however, certain hazards may arise in extraneous circumstances. Do not open the housing of the balance; it does not contain any parts that can be maintained, repaired or replaced by the user. If you experience problems with your balance, contact your authorized METTLER TOLEDO dealer or service representative.

The balance has been tested for the experiments and intended purposes documented in the appropriate manual. However, this does not absolve you from the responsibility of performing your own tests of the products supplied by us regarding their suitability for the methods and purposes you intend to use them for.

#### Intended use

This balance is designed to be used in analytical laboratories by qualified staff. Your balance is used for weighing. Use the balance exclusively for this purpose.

Any other type of use and operation beyond the limits of technical specifications without written consent from Mettler-Toledo GmbH, is considered as not intended.

#### Site requirements

The balance has been developed for indoor operation. Avoid the following environmental influences:

- Conditions outside of the ambient conditions specified in the technical data
- Powerful vibrations
- Direct sunlight
- · Corrosive gas atmosphere

Safety Information Jewelry Balances

- Explosive atmosphere of gases, steam, fog, dust and flammable dust
- Powerful electric or magnetic fields

#### Staff qualification

Incorrect use of the balance or the chemicals used in the analysis can lead to death or injury. The following experience is needed for operating the balance.

- Knowledge and experience in working with toxic and caustic substances.
- Knowledge and experience in working with standard laboratory equipment.
- Knowledge and experience in working in accordance with general lab safety rules.

#### Responsibilities of the balance owner

The balance owner is the person that uses the balance for commercial use or places the balance at the disposal of his staff. The balance owner is responsible for product safety and the safety of staff, user(s) and third party.

The operator has the following responsibilities:

- Know the rules for safety at the workplace that are in effect and enforce them.
- Ensure that only qualified staff uses the balance.
- Define the responsibilities for installation, operation, cleaning, troubleshooting and maintenance and ensure that the tasks are done.
- Train the staff in regular intervals and inform them about dangers.
- · Provide the necessary protective gear for the staff.

#### Shut down of the balance in emergency situations

- Pull the plug from the electrical outlet.

#### Safety notes



### **MARNING**

### Danger of death or serious injury due to electric shock!

Contact with parts that contain a live current can lead to injury and death. If the balance cannot be shut down in an emergency situation, people can be injured or the balance can be damaged.

- 1 Only use the supplied three-core power cable with equipment grounding conductor to connect your balance.
- 2 Check that the voltage printed on the balance is the same as your local power supply voltage.
  - ⇒ If this is not the case, under no circumstances connect the AC/DC adapter to the power supply, but contact a METTLER TOLEDO representative.
- 3 Only connect the balance to a three-pin power socket with earthing contact.
- 4 Only standardized extension cable with equipment grounding conductor must be used for operation of the balance.
- 5 Do not disconnection the equipment grounding conductor.
- 6 Check the cables and the plug for damage and replace damaged cables and plugs.
- 7 Make sure that the cables are arranged so that they cannot be damaged or interfere with the operation.

5

- 8 Keep all electrical cables and connections away from liquids.
- 9 Make sure that the power plug is accessible at all times.



### **NOTICE**

#### **Environment**

Only use indoors in dry locations.

Jewelry Balances Safety Information



# **NOTICE**

### Danger of damaging the keyboard with pointed or sharp objects!

Do not use pointed or sharp objects to navigate on the keyboard. This may damage the surface of the keyboard.

Operate the keyboard with your fingers.



# **NOTICE**

#### Danger of damage to the balance!

Never open the balance. The balance contains no user-serviceable parts.

- In the event of problems, please contact a METTLER TOLEDO representative.



# **NOTICE**

### Danger of damage to the balance due to incorrect parts!

Using incorrect parts with the balance can damage the balance or cause the balance to malfunction.

 Only use parts supplied with the balance, listed accessories and spare parts from Mettler-Toledo GmbH.



# NOTICE

### Damage of the balance or software

In some countries, excessive mains voltage fluctuations and strong glitches may occur. This may affect the balance functions or damage the software.

- Use a voltage regulator for stabilizing.

#### Finding more information




Refers to an external document.

Safety Information Jewelry Balances

# 3 Design and Function

# 3.1 Overview

# 3.1.1 Components

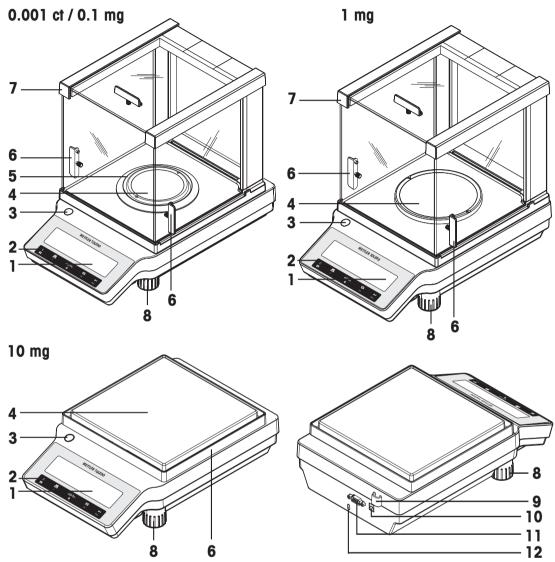


1	Display	2	Operation keys
3	Level indicator	4	Weighing pan
5	Draft shield element	6	Handle for operation of the draft shield door
7	Glass draft shield	8	Leveling foot
9	Legal for Trade (LFT) sealing	10	Socket for AC adapter
11	RS232C serial interface	12	Kensington slot for anti-theft purposes

7

Jewelry Balances Design and Function

# 3.1.2 Operation keys

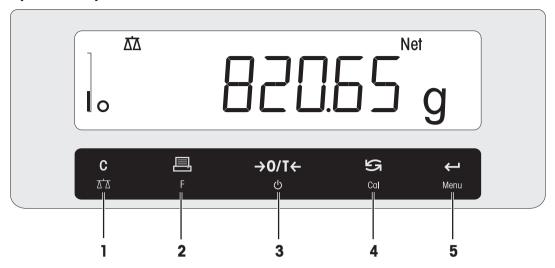


# Legend key functions

8

No.	Key	Press briefly (less than 1.5 s)	Press and hold (longer than 1.5 s)
1	C	<ul><li>Cancel or leave menu without saving</li><li>One step back in the menu</li></ul>	<ul><li>Select the simple weighing application</li><li>Exit application</li></ul>
2	F	<ul> <li>Print display value</li> <li>Transmit data</li> <li>To navigate backwards in the menu or menu selection</li> <li>Decrease parameters in menu or appli- cations</li> </ul>	Open the application list for selecting an application
3	<b>→</b> 0/T← ტ	<ul><li>Zero/Tare</li><li>Switch on</li></ul>	Switch off into standby mode
4	<b>∑</b> Cal	<ul> <li>With entries, scroll down</li> <li>To navigate forward menu topics or menu selections</li> <li>To toggle between unit 1, recall value (if selected), unit 2 (if different from unit 1) and the application unit (if any)</li> <li>Increase parameters in menu or applications.</li> </ul>	Select adjustment (calibration)  with internal weight *  with external weight  Customer fine adjustment *  * On models with internal weight only
5	<b>—</b> Menu	<ul> <li>Enter or leave menu selection</li> <li>To enter application parameter digit and switch to next parameter digit</li> <li>To accept parameter in menu selection.</li> </ul>	<ul> <li>Enter or leave menu (parameter settings)</li> <li>To store parameter</li> <li>To accept numeric inputs in applications.</li> </ul>

Design and Function Jewelry Balances

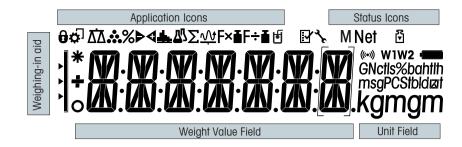

Applica	Application Icons						
$\Delta \Delta$	Application weighing	Σ	Application totaling				
***	Application piece counting	F×∎	Application multiplication factor				
%	Application percent weighing	F÷∎	Application division factor				
<b>▶</b> ∢	Application check weighing	þ	Application density				
<u>.II</u>	Application statistics	0	Menu locked				

#### Note

While an application is running, the corresponding application icon appears at the top of the display.

Status Icons							
M	Indicates stored value (Memory)	(((•))	Acoustic feedback for pressed keys activated				
¥	Adjustments (calibration) started	W1	Weighing range 1 (Dual Range models only)				
3.	Service reminder	W2	Weighing range 2 (Dual Range models only)				

Weight	Weight Value Field and Weighing-in aid							
_	Indicates negative values		Brackets to indicate uncertified digits (approved models only)					
0	Indicates unstable values		Marking of nominal or target weight					
*	Indicates calculated values		Marking of tolerance limit T+					
		Þ	Marking of tolerance limit T-					

Unit Field								
GNctls%bahtlh	g	gram	ozt	troy ounce	tls	Singapore taels		
msgPCStbldigit	kg	kilogram	GN	grain	tit	Taiwan taels		
kgmgm	mg	milligram	dwt	pennyweight	tola	tola		
	ct	carat	mom	momme	baht	baht		
	lb	pound	msg	mesghal				
	OZ	ounce	tlh	Hong Kong taels				


9

Jewelry Balances Design and Function

# 3.2 Basic principles for operation

#### Selecting simple weighing or terminate application

- Press and hold A until WEIGH appears on the display.
  - ⇒ The balance returns to the simple weighing mode.

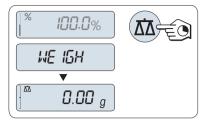


#### Note

How to perform simple weighing **see** [Performing a simple weighing ▶ Page 22].

#### Selecting an application

- Press and hold F until APP.LIST (application list).
   ⇒ Last active application e.g. COUNT appears on the display.
- 2 Select an application by multiple pressing 5.
- 3 To execute selected application press  $\leftarrow$   $\bot$ .

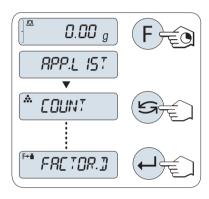


### **Available applications**

Display	Remark	Description
COUNT	Piece counting	see [Application piece counting ▶ Page 34]
PERCENT	Percent weighing	see [Application percent weighing ▶ Page 37]
CHECK	Check weighing	see [Application check weighing ▶ Page 39]
STAT	Statistics	see [Application statistics ▶ Page 41]
TOTAL	Totaling	see [Application totaling ▶ Page 43]
FACTOR.M	Multiplication factor	see [Application multiplication factor weighing ▶ Page 44]
FACTOR.D	Division factor	see [Application division factor weighing ▶ Page 46]
DENSITY	Density	see [Application density ▶ Page 48]

### **Entering the menu**

- 1 Press and hold **Menu** to enter main menu. The first menu **BASIC** is displayed (except menu protection is active).
- 2 Press 🗲 repeatedly to change menu.
- 3 Press ← to confirm the selection.

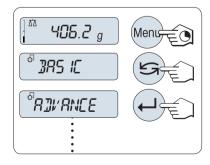

#### Note


10

Detailed description of the menu **see** [The Menu ▶ Page 25].

Design and Function Jewelry Balances


# Selecting menu topics

Press . The next menu topic appears in the display. Each time
 is pressed, the balance switches to the next menu topic.

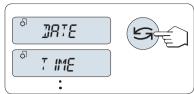

#### Changing settings in selected menu topic

- 1 Press —I. The display shows the current setting in the selected menu topic. Each time is pressed, the balance switches to the next selection. After the last selection, the first is shown again.

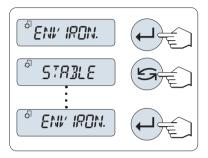


#### Changing settings in a submenu selection

The same procedure as for menu topics.


#### Input principle of numerical values

- Press I to select a digit (cyclically from left to right) or a value (depending on the application). The selected digit or the selected value is blinking.
- 2 For changing blinking digits or values, press at to increase or **F** to decrease.
- 3 Press and hold \leftarrow to accept the value.

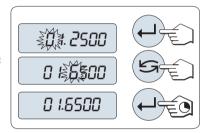

#### Saving settings and closing the menu

- 1 Press and hold **Menu** to leave menu topic.
  - ⇒ **SAVE:YES** appears on the display.
- 2 Press of to toggle between **SAVE:YES** and **SAVE:NO**.
- 3 Press to execute **SAVE:YES**. Changes are saved.
- 4 Press to execute **SAVE:NO**. Changes are not saved.

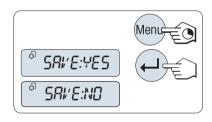

#### Cancel

- During menu operation
- To leave menu topic or menu selection without saving press C (one step back in the menu).
- During application operation
- To cancel settings press C.
  - ⇒ The balance returns to the previous active application.



11

#### **Important**

If no entry is made within 30 seconds, the balance reverts to last active application mode. Changes are not saved. If changes are made, the balance asks "SAVE:NO".

Jewelry Balances Design and Function

# 4 Installation and Putting into Operation



# **WARNING**

### Danger of death or serious injury due to electric shock!

The balance must be disconnected from the power supply when carrying out all setup and mounting work.

# 4.1 Unpacking and delivery inspection



# **NOTICE**

### Danger of damage to the balance due to incorrect parts!

Using incorrect parts with the balance can damage the balance or cause the balance to malfunction.

- Only use parts supplied with the balance, listed accessories and spare parts from Mettler-Toledo GmbH.
- 1 Open the packaging and carefully remove all components.
- 2 Check the delivered items.

#### The standard scope of delivery contains the following items:

Components		Carat balances	Gold balances	Gold balances	
		0.001 ct / 0.1 mg	1 mg	10 mg	
Draft shield	170 mm	✓	✓	_	
Weighing pan	ø 90 mm	✓	_	_	
	ø 120 mm	_	<b>√</b>	_	
	180 × 180 mm	_	_	✓	
Draft shield element		✓	_	✓	
Pan support		_	_	✓	
Protective cover		✓	✓	<b>√</b>	
Universal AC adap	oter (with plug set)	✓	✓	✓	
Carat pan	S 80 ø × 20 mm	✓	_	_	
	M 90 ø × 30 mm	_	✓	_	
EC declaration of conformity		✓	✓	<b>√</b>	
Operating instructions or User Manual; printed or on CD-ROM, depending on country of use		<b>✓</b>	<b>✓</b>	/	

# 4.2 Installing components


## Carat balances with readability of 0.001 ct / 0.1 mg

Place the following components on the balance in the specified order:

- 1 Push the side glass doors back as far as will go.
- 2 Place draft shield element (1).
- 3 Place weighing pan (2).

#### Note

Cleaning the draft shield **see** [Cleaning and service ▶ Page 57].

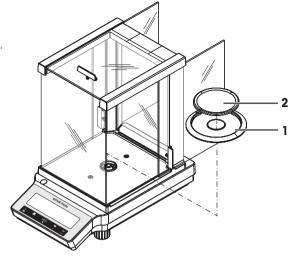


## Gold balances with readability of 1 mg

Place the following components on the balance in the specified order:

- 1 Push the side glass doors back as far as will go.
- 2 Place weighing pan (1).

#### Note


Cleaning the draft shield **see** [Cleaning and service ▶ Page 57].

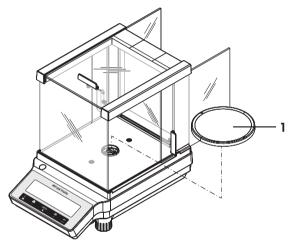

# Gold balances with readability of 10 mg

Place the following components on the balance in the specified order:


- 1 Place draft shield element (1): carefully pull apart the draft shield element to fix it under the retaining plate.
- 2 Insert pan support (2).
- 3 Place weighing pan (3).

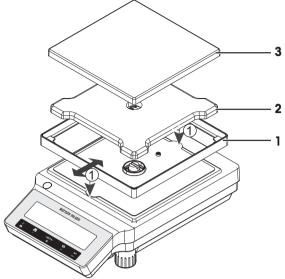

# 4.3 Installing protective cover

### **Important**

Make sure using the correct protective cover, **see** [Accessories and Spare Parts ▶ Page 72].

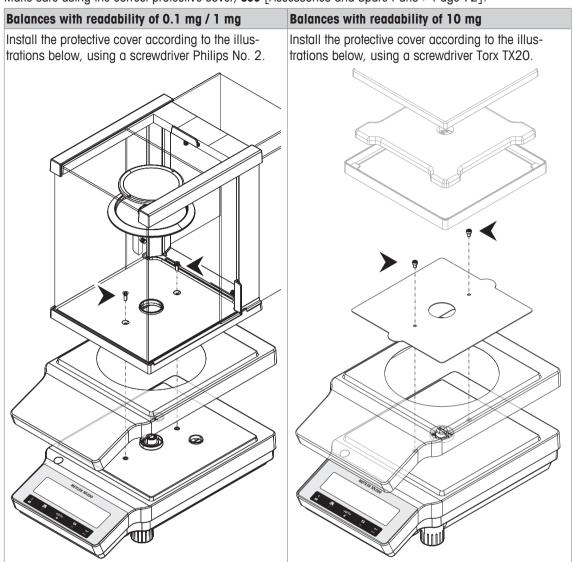

# 4.4 Selecting a location

An optimal location will ensure accurate and reliable operation of the balance. The surface must be able to safely take the weight of the balance when fully loaded. The following local conditions must be observed:

#### **Important**

If the balance is not horizontal at the outset, it must be leveled during commissioning.

- The balance must only be used indoors and up to a maximum altitude of 4,000 m above sea level.
- Before switching on the balance, wait until all parts are at room temperature (+5 to 40 °C).
  - The humidity must be between 10% and 80% non-condensing.
- The power plug must be accessible at all times.
- Firm, horizontal and vibration-free location.
- Avoid direct sunlight.
- No excessive temperature fluctuations.
- No strong drafts.









# 4.5 Connecting the balance



# **↑** WARNING

# Danger of death or serious injury due to electric shock!

Contact with parts that contain a live current can lead to injury and death. If the balance cannot be shut down in an emergency situation, people can be injured or the balance can be damaged.

- 1 Only use the supplied three-core power cable with equipment grounding conductor to connect your balance.
- 2 Check that the voltage printed on the balance is the same as your local power supply voltage.
  - ⇒ If this is not the case, under no circumstances connect the AC/DC adapter to the power supply, but contact a METTLER TOLEDO representative.
- 3 Only connect the balance to a three-pin power socket with earthing contact.
- 4 Only standardized extension cable with equipment grounding conductor must be used for operation of the balance.
- 5 Do not disconnection the equipment grounding conductor.
- 6 Check the cables and the plug for damage and replace damaged cables and plugs.
- 7 Make sure that the cables are arranged so that they cannot be damaged or interfere with the operation.
- 8 Keep all electrical cables and connections away from liquids.
- 9 Make sure that the power plug is accessible at all times.



# **NOTICE**

### Danger of damage to the AC/DC adapter due to overheating!

If the AC/DC adapter is covered or in a container, it is not sufficiently cooled and overheats.


- 1 Do not cover the AC/DC adapter.
- 2 Do not put the AC/DC adapter in a container.

The balance is supplied with an universal AC adapter and a country-specific plug. The AC adapter is suitable for use with the following voltage range:

100 - 240 V AC, 50/60 Hz.

- Connect the AC adapter to the connection socket on the back of your balance (see figure) and to the power line.
  - The balance performs a display test (all segments in the display light up briefly), WELCOME, Software version, Maximum load and Readability appears briefly.

The balance is ready for use.

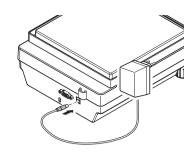

# 4.6 Setting up the balance

### 4.6.1 Switching on the balance

Before working with the balance, it must be warmed up in order to obtain accurate weighing results. To reach operating temperature, the balance must be connected to the power supply for at least 30 minutes (0.1 mg models 60 minutes).

#### Switching on

- The balance is in STANDBY mode. MT.GREEN appears on the display.
- Press (b) or remove any load from weighing pan or tap on the weighing pan.

The balance is ready for weighing or for operation with the last active application.



Approved balances can only be switched on by pressing  $\circlearrowleft$  in selected countries.



# 4.6.2 Leveling the balance

#### **Important**

If the balance is not horizontal at the outset, it must be leveled during commissioning.


### **Important**

The balance must be leveled and adjusted each time it is moved to a new location.

- 1 Align the balance horizontally.
- 2 Turning the two front leveling screws of the housing until the air bubble is in the inner circle of the level indicator.
  - ⇒ The position of the air bubble illustrates which leveling screw you need to turn (L = left leveling screw, R = right leveling screw) and in which direction so that the air bubble moves to the center.



In this example, turn the left leveling screw counterclockwise.

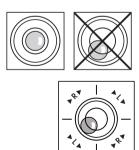


# 4.6.3 Setting date and time

When you put your new instrument into operation for the first time, you should enter the current date and time.

#### Note

- These settings are retained even if you disconnect your instrument from the power supply.
- A reset of the instrument will not change these settings.

- Set the current date according to the date format DATE.FRM in the menu ADVANCE..
- Set the current time according to the time format TIME.FRM in the menu ADVANCE...
- 1 Press and hold Menu until menu BASIC appears on the display.
- 2 Press ← to open menu BASIC.
  - ⇒ DATE appears.
- 3 Press ← to confirm.
- 4 **Set current date**. Press ← to select day, month or year; press to set current day, month or year.
- 5 Press and hold ← to confirm the settings.
  - ⇒ DATE appears...
- 6 Set current time. Press 🗲 to select TIME.
- 7 Press to confirm.
  - ⇒ +1H appears.
- 8 Select **SET.TIME** by pressing **S**.
- 9 Press ← to confirm.
- 11 Press and hold \(\rightarrow\) to confirm the settings.
  - ⇒ **TIME** appears.
- 12 Press and hold \to store the settings.
  - ⇒ **SAVE:YES** appears.
- 13 Press ← to confirm.

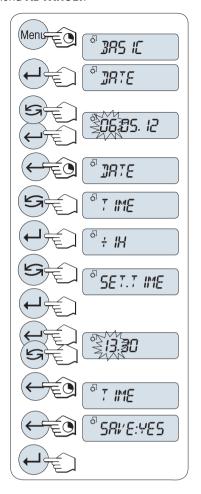

# 4.6.4 Adjusting the balance

To obtain accurate weighing results, the balance must be adjusted to match the gravitational acceleration at its location and depending on the ambient conditions. After reaching the operation temperature, adjusting is necessary

- before the balance is used for the first time.
- after a change of the location.
- at regular intervals during weighing service.

# 4.7 Adjustment (calibration)



### NOTICE

Before adjusting the balance, it must be warmed up.


# 4.7.1 Adjustment with internal weight

#### Note

On models with internal weight only (see technical data).

- Weighing pan is unloaded.
- To carry out this operation press and hold CAL until ADJUST appears.
- 2 Select **ADJ.INT** by pressing **S**.
  - ⇒ **ADJ.INT** appears on the display.
- 3 Press ← to execute internal adjustment.

The balance adjusts itself automatically. The adjusting is finished when the message **ADJ.DONE** appears briefly on the display. The balance returns to the last active application and is ready for operation.

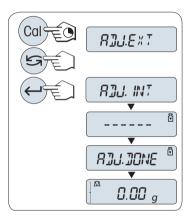

### Sample adjustment printout using internal weight:

```
- Internal Adjustment --
21.Jan 2012 12:56

METTLER TOLEDO

Balance Type JE703C
SNR 1234567890

Temperature 22.5 °C
Diff 3 ppm


Adjustment done
```

# 4.7.2 Adjustment with external weight

#### **Important**

Because of certification legislation, the approved models cannot be adjusted with an external weight \* (depend on selected countries' certification legislation).

- \* except OIML accuracy class I approved models.
- 1 Have required adjustment weight ready.
- 2 To carry out this operation press and hold CAL until ADJUST appears.
- 3 Select **ADJ.EXT** by pressing **S**.
  - ⇒ **ADJ.EXT** appears on the display.
- 4 Unload weighing pan.
- 5 Optional: If necessary, you can define a different weight value. Press to change a digit (cyclically from left to right); press to change the blinking digit.
- 6 Press and hold \(\rightarrow\) to execute external adjustment.
  - ⇒ The required adjustment weight value flashes in the display.
- 7 Place adjustment weight in center of pan.
  - ⇒ The balance adjusts itself automatically.
- 8 When zero is flashing, remove adjustment weight.
- ⇒ The adjusting is finished when the message ADJ.DONE appears briefly on the display. The balance returns to the last active application and is ready for operation.

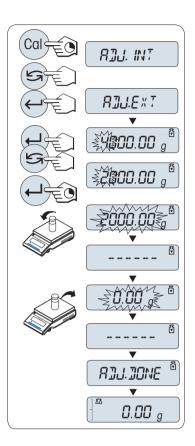

#### Sample adjustment printout using external weight:

```
- External Adjustment --
21.Jan 2012 12:56

METTLER TOLEDO

Balance Type JE3002GE
SNR 1234567890

Temperature 22.5 °C
Nominal 2000.00 g
Actual 1999.99 g
Diff 5 ppm

Adjustment done

Signature
```

### 4.7.3 Customer fine adjustment



# **NOTICE**

This function should be executed only by trained personnel.

The function customer fine adjustment **ADJ.CF** allows you to adjust the value of the internal adjustment weight with your own adjustment weight. The adjustable range of the adjustment weight is possible only in a very small range. Customer fine adjustment impacts the function of internal adjustment. The customer fine adjustment can be deactivated at any time.

#### Note


- This feature is available on models with internal weight only.
- Because of certification legislation, approved models cannot be adjusted with customer fine adjustment (depending on selected countries' certification legislation).
- Use certificated weights.
- Balance and test weight have to be on operating temperature.
- Observe the correct environmental conditions.

#### **Execute customer fine adjustment**

- The balance is under measuring condition.
- 1 Have required adjustment weight ready.
- 2 Unload weighing pan.
- 3 To carry out this operation press and hold CAL until ADJUST appears.
- 4 Select **ADJ.CF** by pressing **S**.
  - ⇒ ADJ.CF appears on the display.
- 5 Select **EXECUTE**.
- 6 Start Adjustment with ← J.
  - ⇒ **SET REF.** appears briefly.
  - ⇒ The last saved value flashes on the display.
- 7 Select the target adjustment weight. Press ← to change a digit (cyclically from left to right); press ← to change the blinking digit.
- 8 Press and hold \(\llorer\) to confirm and execute **ADJ.CF**.
  - ⇒ The required adjustment weight value flashes in the display. This could take some time.
- 9 Place required adjustment weight in center of pan.
- 10 Remove adjustment weight when zero is flashing.
- 11 Wait until ADJ.DONE briefly appears.
- ⇒ The adjusting is finished when the message ADJ.DONE appears briefly on the display. The balance returns to the last active application and is ready for operation.
- ⇒ If the error message WRONG ADJUSTMENT WEIGHT appears, the weight is not within the allowed value range and could not be accepted. ADJ.CF could not be executed.

#### Note

Storing the adjustment is not required.

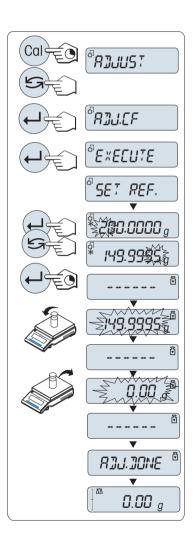


#### Deactivate customer fine adjustment

- 1 To carry out this operation press and hold **CAL** until **ADJUST** appears.
- 2 Select **ADJ.CF** by pressing **S**.
  - ⇒ **ADJ.CF** appears on the display.
- 3 Select RESET.
- 4 Start **RESET** by pressing ← ...
  - ⇒ NO? appears.
- 5 Select YES? and confirm with ← .
- ⇒ The adjusting is finished when the message **ADJ.DONE** appears briefly on the display. The balance returns to the last active application and is ready for operation with initial adjustment.

# 4.8 Performing a simple weighing


- 1 Press  $\rightarrow 0/T \leftarrow$  to zero the balance.
  - **Note:** If your balance is not in the weighing mode, press and hold the  $\overline{\Delta\Delta}$  key down until **WEIGHING** appears in the display. Release the key. Your balance is in the weighing mode and set to zero.
- 2 Place weighing sample on the weighing pan.
- 3 Wait until the instability detector o disappears and the stability beep sounds.
- 4 Read the result.

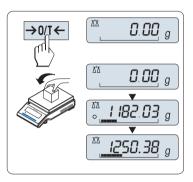

#### Zeroing


- 1 Unload the balance.
- 2 Press → 0/T ← to set the balance to zero. All weight values are measured in relation to this zero point.

Use the  $\rightarrow 0/T \leftarrow$  zeroing key before you start with a weighing.

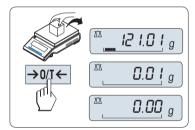


#### **Taring**

- If you are working with a weighing container, first set the balance to zero
- Place empty container on the weighing pan. The weight is displayed.
- 2 Press  $\rightarrow 0/T \leftarrow$  to set the balance to zero.
  - ⇒ **0.00 g** appears in the display.
- 3 Place weighing sample into the weighing container.
- ⇒ The result appears in the display.

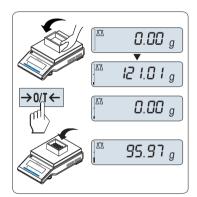

#### Switching weight units

The key can be used at any time to toggle between weight unit **UNIT 1**, **RECALL** value (if selected) and weight unit **UNIT 2** (if different from weight unit 1) and the application unit (if any).

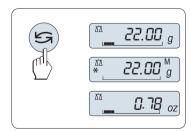

#### Recall / Recall weight value

Recall stores stable weights with an absolute display value bigger than 10d.

**Requirement:** The function **RECALL** must be activated in the menu.

- 1 Load weighing sample. The display shows weight value and stores stable value.
- 2 Remove weighing sample. When the weight is removed the display shows zero.
- 3 Press . The display shows last stored stable weight value for 5 seconds together with asterisk (\*) and memory (M) symbols. After 5 seconds the display goes back to zero. This can be repeated unlimited times.

#### Delete last weight value

As soon a new stable weight value is displayed, the old recall value becomes replaced by the new weight value. When pressing  $\rightarrow 0/T \leftarrow$ , the recall value is set to 0.

**Note:** If the power is switched off, the recall value is lost. The recall value can not be printed.

#### Weighing with the weighing-in aid

The weighing-in aid is a dynamic graphic indicator which shows the used amount of the total weighing range. You can thus recognize at a glance when the load on the balance approaches the maximum load.

# 0% 100%

0.00 a

12 1.0 1 g

0.00 a

 $0.00_{g}$ 

12 1.0 1

#### Print / Transmit data

Pressing the 🗏 key transmits the weighing results over the interface e.g. to a printer or a PC.

### Switching off into standby mode

- Press and hold the key until **STANDBY** appears on the display.
   Release the key.
- ⇒ MT.GREEN appears on the display.





#### Note

- Once your balance has been switched off, it is in energy saver mode **STANDBY**. In this case your balance needs no warm-up time in the standby mode and is immediately ready for weighing. If you wish to perform a weighing, you only need to place the sample on the weighing pan and the balance immediately displays the result. There is no need to switch it on with the **b** key (with approved balances only possible in selected countries).
- To completely switch off the balance, disconnect it from the power supply.

# 4.9 Weighing below the balance

Your balance is equipped with a hanger for carrying out weighings below the work surface (weighing below the balance).




# **NOTICE**

#### Damage to balance

Do not place the balance on the pan support location bolt.

- 1 Press and hold the  $\bigcirc$  key.
- 2 Disconnect the balance from the power supply.
- 3 Disconnect all interface cables.
- 1 Remove weighing pan, pan support and draft shield element if present.
- 2 Turn the balance carefully on its side.
- 3 Remove the cap. Keep it for later use.
- 4 Turn the balance to its normal position and simply reinstall all components in the reverse order.

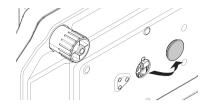

# 4.10 Transporting the balance

### Switching off the balance

- 1 Press and hold the 🖒 key.
- 2 Disconnect the balance from the power supply.
- 3 Disconnect all interface cables.

#### Transport over short distances

To move the balance over a short distance to a new location, follow the instructions below.



### NOTICE

### Danger of damage to the glass draft shield!

Glass parts of the balance can be damaged.

 Never lift the balance by the glass draft shield. The draft shield is not sufficiently fastened to the balance.

#### Transporting over long distances

The complete original packaging must be used for transportation or shipment of the balance over long distances or if it cannot be ensured that the balance will be transported upright.

# 5 The Menu

# 5.1 What is in the menu?

The menu allows you to match your balance to your specific weighing needs. In the menu you can change the settings of your balance and activate functions. The main menu has 4 different menus and these contains different topics, each of which allows you various selection possibilities.

For menu **PROTECT see** [Main menu ▶ Page 26].

#### **Menu BASIC**

Topic	Explanation	Description
DATE	Setting the current date.	[see ▶ Page 26]
TIME	Setting the current time.	[see ▶ Page 26]
1/10 D	Setting display increment (1/10d function)	[see ▶ Page 26]
UNIT 1	Specification of the 1 <sup>st</sup> weight unit in which the balance should show the result.	[see ▶ Page 26]
UNIT 2	Specification of the 2 <sup>nd</sup> weight unit in which the balance should show the result.	[see ▶ Page 27]
SET ID	Setting an identification.	[see ▶ Page 27]
PRT.MENU	Printing the settings.	[see ▶ Page 27]
RESET	Call up of the factory settings.	[see ▶ Page 27]

#### Menu ADVANCE.

Topic	Explanation	Description
ENVIRON.	Matching the balance to the ambient conditions.	[see ▶ Page 27]
ADJ.LOCK	Switching the adjustment function on or off.	[see ▶ Page 28]
DATE.FRM	Setting the date format.	[see ▶ Page 28]
TIME.FRM	Preselection of the time format.	[see ▶ Page 28]
RECALL	Switching the application "Recall" for storing stable weights on or off.	[see ▶ Page 28]
STANDBY	Setting the time after which the balance should be switched off automatically.	[see ▶ Page 28]
B.LIGHT	Switching on or off the display backlight.	[see ▶ Page 28]
A.ZERO	Switching the automatic zero correction (Autozero) on or off.	[see ▶ Page 29]
SRV.ICON	Switching the service reminder (service icon) on or off.	[see ▶ Page 29]
SRV.D.RST	Reset service date and hours (service reminder)	[see ▶ Page 29]

#### **Menu INT.FACE**

Topic	Explanation	Description
RS232	Matching the serial interface RS232C to a peripheral unit.	[see ▶ Page 29]
HEADER	Setting the header for printout of individual values.	[see ▶ Page 30]
SIGN.L	Setting the footer for printout of individual values.	[see ▶ Page 30]
LN.FEED	Setting line feed for printout of individual values.	[see ▶ Page 30]
ZERO.PRT	Setting the auto print function for printing zero.	[see ▶ Page 30]
COM.SET	Setting the data communication format of the serial interface RS232C.	[see ▶ Page 30]
BAUD	Setting the transfer speed of the serial interface RS232C.	[see ▶ Page 31]
BIT.PAR.	Setting the character format (Bit/Parity) of the serial interface RS232C.	[see ▶ Page 31]

25

Jewelry Balances The Menu

Topic	Explanation	Description
STOPBIT	Setting the character format (stop bit) of the serial interface RS232C.	[see ▶ Page 32]
HD.SHK	Setting the transfer protocol (Handshake) of the serial interface RS232C.	[see ▶ Page 32]
RS.TX.E.O.L.	Setting the end of line format of the serial interface RS232C.	[see ▶ Page 32]
RS.CHAR	Setting the char set of the serial interface RS232C.	[see ▶ Page 32]
INTERVL.	Selection of the time interval for the simulated print key press.	[see ▶ Page 32]

# 5.2 Description of menu topics

In this section you will find information regarding the individual menu topics and the available selections.

### 5.2.1 Main menu

Selecting the submenu.

**BASIC** The small **BASIC** menu for simple weighing is displayed. **ADVANCE.** The extended **ADVANCE.** menu for further weighing settings is

displayed.

**INT.FACE** The menu **INT.FACE** for all interface parameter settings for

peripheral devices e.g. printer is displayed.

**PROTECT** Menu protection. Protection of balance configurations against

unmeant manipulation.

**OFF** Menu protection is off. (**Factory setting**)

**ON** Menu protection is on. The menu **BASIC**, **ADVANCE**. and

**INT.FACE** are not displayed. This is indicated with **1** in the

display.

#### 5.2.2 Basic menu

#### DATE - Date

Setting the current date according to date format.

**Important:** A reset of the balance will not change this setting.

#### TIME - Time

Setting the current time according to time format

**+1H** Set the current time forwards by 1 hour. (Factory setting)

**-1H** Set the current time backwards by 1 hour.

**SET.TIME** Enter the current time.

**Important:** A reset of the balance will not change this setting.

#### 1/10 D - Display increment 1/10 d

This menu topic allows you to reduce the readability of the display.

**Important:** This menu topic is not available with models which are approved and e=d.

**OFF** 1/10 **D** display increment is switched off (full resolution)

(Factory setting)

**ON** 1/10 **D** switched on (low resolution)

Important: A reset of the balance will not change this setting.

#### UNIT 1 - Weight unit 1

Depending on requirements, the balance can operate with the following units (depending on the model)

26 | The Menu Jewelry Balances

- Only those weight units allowed by the appropriate national legislation are selectable.
- With approved balances, this menu topic has a fixed setting and cannot be changed.

g	Gram	dwt	Pennyweight
kg	Kilogram	mom	Momme
mg	Milligram	msg	Mesghal
ct	Carat	tlh	Tael Hong Kong
lb	Pound	tis	Tael Singapore
OZ	Ounce (avdp)	tlt	Tael Taiwan

tola

baht

#### UNIT 2 - Weight unit 2

Units:

ozt

GN

If it is required to show the weighing results in weighing mode in an additional unit, the desired second weight unit can be selected in this menu topic (depending on the model). Units see UNIT 1.

**Important:** Only those weight units allowed by the appropriate national legislation are selectable.

#### SET ID - Set identification

This menu topic allows you to set your own desired identification to the balance for the convenience of asset management or other purposes. The ID can be printed with other balance information. One ID can be set and max 7 alphanumeric characters are possible (blank, 0...9, A...Z).

> **SET ID** Set identification

Ounce (troy)

Grain

The setting starts from left to right and the display prompts the configurable position by flashing corresponding place.

Tola

Baht

- SET ID is selected.
- 1 Search through (blank, 0...9, A...Z) by pressing 5.
- 2 After selecting the character, press ← to confirm and move to the next place. To store press and hold  $\leftarrow$ .

**Important:** A reset of the balance will not change this setting.

#### PRT.MENU - Print menu

This menu topic allows you to execute a printout of the menu settings if a printer is connected. This topic is only visible if **PRINTER** mode is selected.

- PRT.MENU appears on the display and a printer is properly connected.
- To execute a printout press ←

#### **RESET – Reset balance settings**

This menu topic allows you to call-up the factory settings.

To toggle between **YES**? and **NO**? press **S**.

Important: A reset of the balance will not change DATE, TIME, 1/10 D, and SET ID settings.

#### 5.2.3 Advanced menu

#### ENVIRON. – Environment settings

This setting can be used to match your balance to the ambient conditions.

STD. Setting for an average working environment subject to moderate

variations in the ambient conditions. (Factory setting)

27

UNSTAB. Setting for a working environment where the conditions are

continuously changing.

**STABLE** Setting for a working environment which is practically free from

drafts and vibrations.

Jewelry Balances The Menu

#### ADJ.LOCK - Adjustment (calibration) lock

Under this menu topic you can lock function of the Cal key.

**OFF** The adjustment lock is switched **off**. The adjustment function is

on. The Cal key is activ. (Factory setting)

**ON** The adjustment lock is switched **on**. The adjustment function is

off. The Cal key has no function.

#### DATE.FRM - Date format

This menu topic allows you to preselect the date format.

The following date formats are available:

	Display examples	Printing examples
DD.MM.Y	01.02.09	01.02.2009
MM/DD/Y	02/01/09	02/01/2009
Y-MM-DD	09-02-01	2009-02-01
D.MMM Y	1.FEB.09	1.FEB 2009
MMM D Y	FEB.1.09	FEB 1 2009

### Factory setting: DD.MM.Y

#### TIME.FRM - Time format

This menu topic allows you to preselect the time format.

The following date formats are available:

24:MM 12:MM 24.MM 12.MM

3.04 PM

Factory setting: 24:MM

#### RECALL - Recall

This menu topic allows you to switch the **RECALL** function on or off. When it is switched on recall stores the last stable weight if the absolute display value was bigger than 10d.

OFF	RECALL switched of	off. (Factor	y setting)
-----	--------------------	--------------	------------

ON RECALL switched on.

**Important:** The recall value is displayed with an asterisk and cannot be printed.

#### STANDBY – Automatic standby

If the automatic standby function is activated, the balance automatically switches itself after a pre selected time of inactivity into the energy saver mode **STANDBY** (e.g. with no key being pressed and no changes of weight occurring).

**A.OFF** Automatic standby deactivated.

**A.ON** Automatic standby activated (**Factory setting**).

Time in minutes of inactivity for activating standby function.

Setting range: 2...720 minutes.

#### **B.LIGHT - Backlight**

Under this menu topic, the display backlight can be switched off or on.

**B.L. ON** Backlight is always **on**. (**Factory setting**)

**B.L. OFF** Backlight is always **off**.

The Menu Jewelry Balances

#### A.ZERO – Automatic zero setting

This menu topic allows you to switch the automatic zero setting on or off.

**ON A.ZERO switched on** (factory setting). The automatic zero

setting continuously corrects possible variations in the zero point that might be caused through small amounts of contami-

nation on the weighing pan.

OFF A.ZERO switched off. The zero point is not automatically

corrected. This setting is advantageous for special applications

(e.g. evaporation measurements).

**Important:** With approved balances, this setting is not available (only available in selected countries).

#### SRV.ICON - Service reminder

This menu topic allows you to switch the service reminder  $\$  on or off.

ON Service reminder \*\switched on. You will be informed to call

service for recalibration. This will be indicated by the flashing

service icon: 🔧. (Factory setting)

**OFF** Service reminder  $\$  **switched off**.

#### SRV.D.RST - Service date reset

This menu topic allows you to reset service date.

Important: This menu topic is only available if SRV.ICON setting ON was selected.

To toggle between **YES**? and **NO**? press **S**.

#### 5.2.4 Interface menu

#### RS232 - RS232C interface

At this menu topic you can select the peripheral device connected to the RS232C interface and specify how the data is transmitted.

**PRINTER** Connection to a printer. (Factory setting)

Note

Only one printer possible.

**(3)** 

Refer to your printer documentation for recommended printer

settings.

**PRT.STAB** If the Rey is pressed, the next stable weight value will be

printed. (Factory setting)

**PRT.AUTO** Every stable weight value will be printed, without pressing the

昌 key.

PRT.ALL If the 🗏 key is pressed, the weight value will be printed

regardless of stability.

**PC-DIR.** Connection to a PC: the balance can send data (as a

Keyboard) to the PC used for PC applications e.g. Excel.

Note

• The balance sends the weight value without the unit to the

29

PC.

Not available on Win7.

**PRT.STAB** If the \( \bullet \) key is pressed, the next stable weight value will be

sent followed by an enter. (Factory setting)

**PRT.AUTO** Every stable weight value will be sent followed by an enter,

without pressing the 🗏 key.

Jewelry Balances The Menu

PRT.ALL If the 🗏 key is pressed, the weight value will be sent followed

by an enter regardless of stability.

**HOST** Connection to a PC, Barcode Reader etc.: the balance can send

data to the PC and receive commands or data from the PC. **Note** The balance sends the complete MT-SICS answer to the PC (see chapter "MT-SICS interface commands and functions".

**SND.OFF** Send mode switched off. (**Factory setting**)

**SND.STB** If the kev is pressed, the next stable weight value will be

sent.

**SND.CONT** All weight value updates will be sent regardless of stability,

without pressing the \bullet key.

**SND.AUTO** Every stable weight value will be sent, without pressing the  $\square$ 

key.

**SND.ALL** If the \( \exists \) key is pressed, the weight value will be sent

regardless of stability.

2.DISP Connection of an optional auxiliary display unit.

**Important:** The transmission parameters cannot be selected.

Settings are automatically set.

#### HEADER – Options for the printout header of individual values

This menu topic allows you to specify the information that is to be printed at the top of the printout for every individual weighing results (after pressing ).

Important: This menu topic is only available if PRINTER setting was selected.

NO The header is not be printed. (Factory setting)

**DAT/TIM** Date and time are printed.

**D/T/BAL** Date, time and balance information (Balance type, SNR,

Balance ID) are printed. **Note:** Balance ID only if set.

#### SIGN.L – Options for the printout footer for signature line of individual values

This menu topic allows you to set a footer for signature at the bottom of the printout for every individual weighing result (after pressing =).

**Important:** This menu topic is only available if **PRINTER** setting was selected.

**OFF** The signature footer is not be printed. (**Factory setting**)

**ON** The signature footer is printed.

#### LN.FEED – Options for complete the printout of individual values

This menu topic allows you to specify the number of blank lines to complete the printout (line feed) for every individual weighing result (after pressing =).

Important: This menu topic is only available if PRINTER setting was selected.

Possible numbers of blank lines: 0 to 99. (Factory setting =

0)

#### **ZERO.PRT – Options for PRT.AUTO**

This menu topic allows you to specify the auto print function PRT.AUTO for printing zero YES or NO.

**OFF** Zero is not be printed (Zero +/- 3d). (**Factory setting**)

**ON** Zero is always printed.

Important: This menu topic is only available if PRT.AUTO function of the PRINTER or PC-DIR. was selected.

#### COM.SET – Options for the data communication format (RS232C) (HOST)

This menu topic allows you to set the data format depending on which peripheral device is connected.

Important: This menu topic is only available if HOST setting was selected.

30 | The Menu Jewelry Balances

MT-SICS The MT-SICS data transfer formats is used. (Factory setting)

For more information see section "MT-SICS interface commands

and functions".

**SART** The following Sartorius commands are supported:

K Ambient conditions: very stable
 L Ambient conditions: stable
 M Ambient conditions: unstable
 N Ambient conditions: very unstable

O Block keys

P Print key (print, auto print; activate or block)

R Unblock keys
S Restart/self-test

T Tare key

W Calibration/adjustment \*)

Z Internal calibration/adjustment \*\*)

fl\_ Function key (CAL)

s3\_ C key

x0\_ Perform internal calibration \*\*)
 x1\_ Print balance/scale model
 x2 Print weighing cell serial number

x3\_ Print software version

#### **Functionality mapping**

HOST settings: Sartorius printer settings:

**SND.OFF** not applicable

 SND.STB
 manually print with stability

 SND.ALL
 manually print without stability

 SND.CONT
 automatically print without stability

 SND.AUTO
 similar applicable to automatically print

when load is changed

31

#### BAUD - Baud rate RS232C

This menu topic allows you to match the data transmission to different serial RS232C receivers. The baud rate (data transfer rate) determines the speed of transmission via the serial interface. For problem-free data transmission the sending and receiving devices must be set at the same value.

The following settings are available:

600 bd, 1200 bd, 2400 bd, 4800 bd, 9600 bd (Factory setting), 19200 and 38400 bd.

#### **Important**

- Not visible for 2nd display.
- Each device has separate settings.

#### BIT.PAR. - Bit/Parity RS232C

At this menu topic you can set the character format for the attached RS232C serial peripheral device.

**8/NO** 8 data bits/no parity (**Factory setting**)

**7/NO** 7 data bits/no parity

Jewelry Balances The Menu

<sup>\*)</sup> May be inaccessible on verified balances/scales

<sup>\*\*)</sup> Only on models with built-in motorized calibration weight

7/MARK	7 data bits/mark parity
7/SPACE	7 data bits/space parity
7/EVEN	7 data bits/even parity
7/ODD	7 data bits/odd parity

#### **Important**

- Not visible for 2nd display.
- Each device has separate settings.

#### STOPBIT - Stop bits RS232C

At this menu topic you can set the stop bits of the transmitted data to different RS232C serial receivers.

1 BIT 1 Stop bit (Factory setting)

2 BITS 2 Stop bits

#### HD.SHK - Handshake RS232C

This menu topic allows you to match the data transmission to different RS232C serial receivers.

XON.XOFF Software handshake (XON/XOFF) (Factory setting)

RTS.CTS Hardware handshake (RTS/CTS)

**OFF** No handshake

#### **Important**

- Not visible for 2nd display.
- Each device has separate settings.

### RS.TX.E.O.L. - End of line RS232C

At this menu topic you can set the end of line character of the outgoing transmitted data to different RS232C serial receivers.

**CR LF** <CR><LF> Carriage Return followed by Line feed (ASCII-Codes

013 + 010) (Factory setting)

**CR** <CR> Carriage Return (ASCII-Code 013)

**LF** <LF> Line feed (ASCII-Code 010)

**DIR.** is selected)

#### **Important**

- Not visible for 2nd display.
- Each device has separate settings.

#### RS.CHAR - Char set RS232C

At this menu topic you can set the character set of the transmitted data to different RS232C serial receivers.

**IBM.DOS** Char set IBM/DOS (**Factory setting**)

ANSI.WIN Char set ANSI/WINDOWS

#### **Important**

- Not visible for 2nd display.
- Each device has separate settings.

#### INTERVL. - Print key simulation

At this menu topic you can activate a simulation of the key. **INTERVL.** simulates a print key press every x seconds.

32 | The Menu Jewelry Balances

Range: 0 to 65535 seconds

0 sec: Disables the print key simulation

Factory setting: 0 sec

Important: The executed action is according to the configuration of the print key, (see interface setting).

Jewelry Balances The Menu 33


# **6** Applications

# 6.1 Application piece counting




The **Piece counting** application allows you to determine the number of pieces put on the weighing pan. All pieces must be of approximately equal weight, since the number of pieces is determined on the basis of average weight.

- 1 Call-up APP.LIST by pressing and holding F.
- 2 Select application **COUNT** by scrolling with **S**.
- 3 Activate function **COUNT** by pressing ← J.

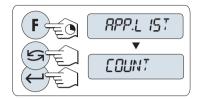


Piece counting first requires the setting of a reference weight, there are 4 possibilities:

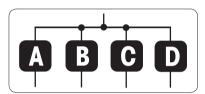

- A Setting the reference by multiple pieces with fix reference values.
- B Setting the reference by multiple pieces with variable reference values.
- C Setting the reference for 1 piece in weighing mode.
- D Setting the reference for 1 piece in manual mode.

Setting the reference by multiple pieces with fix reference values

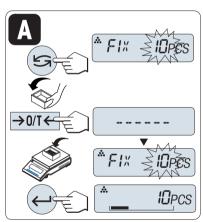

- 1 Select a number of reference pieces by scrolling with S. Possible numbers\* are 5, 10, 20 and 50.


  \* with approved balances in selected countries: min 10.
- 2 Press  $\rightarrow 0/T \leftarrow$  to tare. If using: place empty container on the weighing pan first or tare again.
- 3 Add the selected number of reference pieces to container.
- 4 Press ← to confirm.

Applications Jewelry Balances

#### Setting possibility


# Setting the reference by multiple pieces with variable reference values

- 1 Select **VAR.REF** by scrolling with **⊆**. Press **←** to confirm.
- 2 Select the number of reference pieces. Possible numbers are 1 to 999. With approved balances in selected countries: min 10
- 3 To select a digit, press ← (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 4 To change the digit, press 5.
- 5 Press → **0/T** ← to zero/tare. If using: place empty container on the weighing pan first or zero/tare again.
- 6 Add the selected number of reference pieces to container.
- 7 Press and hold  $\leftarrow$  to confirm.

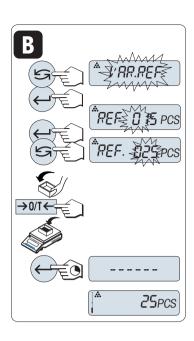

35

#### Setting the reference for one piece in weighing mode

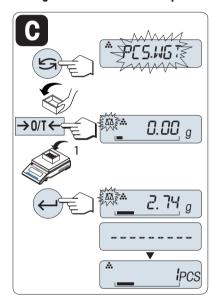

- 1 Select **PCS.WGT** by scrolling with 5.
- 2 Press  $\rightarrow 0/T \leftarrow$  to tare. If using: place empty container on the weighing pan first or tare again.
- 3 Add one reference piece to container. The weight of one piece is displayed.
- 4 Press ← to confirm.

#### **Important**

With approved balances, this setting is not available in selected countries.

Jewelry Balances Applications


#### Setting possibility



#### Setting the reference for one piece in manual mode

- Select **PCS.WGT** by scrolling with **S**.
- 2 Press ← to confirm.
- 3 Enter the final reference one piece weight.
- 4 To select a digit, press ← (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 5 To change the digit, press 5.
- 6 Press and hold ← to confirm.

**Important:** With approved balances, this setting is not available in selected countries.

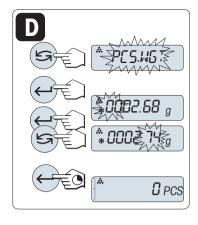

**Important:** If without any key press within 60 seconds, the balance return to the previous active application. Press **C** to cancel and returns to the previous active application.

On completion of the setting procedure, your balance is ready for piece counting.

#### Note

36

- The **RECALL** value is displayed with an asterisk (\*) and icon **M** and can not be printed.
- Take into account minimum values: min. reference weight = 10d (10 digits), min. piece weight\* = 1d (1 digit)!
  - \* with approved balances in selected countries: min 3e
- The current reference weight remains stored until the reference setting is changed.

#### **Exit current application**


To exit the current application and to return to simple weighing mode, press and hold  $\overline{\Delta}\overline{\Delta}$  (longer than 1.5 s).

# 6.2 Application percent weighing




The **Percent weighing** application allows you to check a sample weight as percentage to a reference target weight.

- 1 Call-up **APP.LIST** by pressing and holding **F**.
- 2 Select application **PERCENT** by scrolling with S.
- 3 Activate function **PERCENT** by pressing ← J.

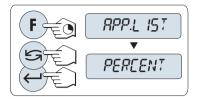

Percent weighing first requires the setting of a reference weight that should corresponds to 100%, there are 2 possibilities:

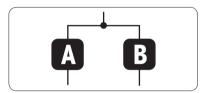


- A Setting the reference in manual mode (enter 100%).
- B Setting the reference in weighing mode (weigh 100%).

Setting possibility




#### Setting the reference by manual mode (enter 100%)

- 1 Press ← to activate manual mode.
- 2 To select a digit, press  $\leftarrow$  (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 3 To change the digit, press 5.
- 4 Press and hold \leftarrow to confirm the value.

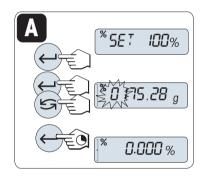

37

#### Setting the reference by weighing mode (weigh 100%)

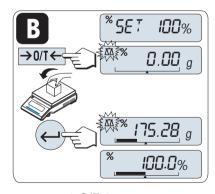


- Press  $\rightarrow 0/T \leftarrow$  to tare the balance and to activate the weighing mode. If needed: place empty container on the weighing pan and tare again.
- 2 Load the reference weight (100%).
  - Note: Reference weight must be at least +/- 10d.
- 3 Press ← to confirm.

**Important:** if without any key press within 60 seconds or by pressing **C**, the balance returns to the previous active application.

On completion of the weighing-in procedure, your balance is ready for percent weighing.

Jewelry Balances Applications

#### Switching between percent and weight display

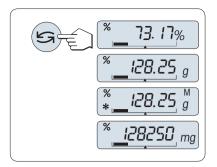

Press key at any time to switch the display between percent display, weighing unit UNIT 1, RECALL value (if activated) and weighing unit UNIT 2 (if different from UNIT 1).

#### Note

38

- The recall value is displayed with an asterisk (\*) as well as icon **M** and can not be printed.
- The current set weight remains stored until it is redetermined.

#### **Exit current application**

To exit the current application and to return to simple weighing mode, press and hold  $\overline{\Delta}\Delta$  (longer than 1.5 s).

# 6.3 Application check weighing



The **Check weighing** application allows you to check the deviation of a sample weight within a tolerance limit to a reference target weight.

- Call-up **APP.LIST** by pressing and holding **F**.
- 2 Select application **CHECK** by scrolling with **S**.
- 3 Activate function **CHECK** by pressing ← .



Step 1: Check weighing first requires the setting of a reference weight that should corresponds to the nominal weight, there are 2 possibilities:

- A Setting the reference in manual mode (enter nominal weight).
- B Setting the reference in weighing mode (weigh nominal weight).

Step 2: Check weighing needs the upper and lower limits:


2 Setting the upper and lower limits in percentage.

Step 1, setting possibility:




- 2 Select the reference target weight.
- 3 To select a digit, press ← (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 4 To change the digit, press 5.
- 5 Press and hold \to confirm the nominal weight.


Step 1, setting possibility:

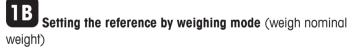


- 1 Press  $\rightarrow 0/T \leftarrow$  to tare the balance and to activate the weighing mode. If using: place empty container on the weighing pan first or tare again.
- 2 Load the nominal weight.
- 3 Press ← to confirm the nominal weight.



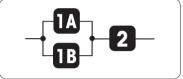

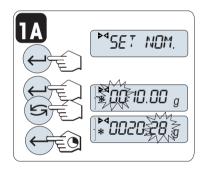

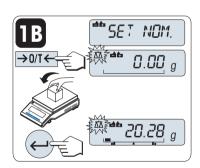

39

**Jewelry Balances Applications** 



# Setting the upper and lower limits (in percentage):

- 1 Press ← to start setting.
- 2 Press ← to confirm the default limit of +/- 2.5% or enter the limit value.
- 3 To select a digit, press  $\longleftarrow$  (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 4 To change the digit, press 5.
- 5 Press and hold ← to confirm the limits.

#### **Important**


- If without any key press within 60 seconds, the balance returns to the previous active application. Press **C** to cancel and returns to the previous active application.
- The nominal weight must be at least 10 digit.

#### On completion of the setting procedure, your balance is ready for checkweighing.

#### Weighing-in-Aid

The Weighing-in-Aid helps you quickly determine the position of the sample weight regarding the tolerance.

- 1 Lower limit
- 2 Target weight
- 3 Upper limit

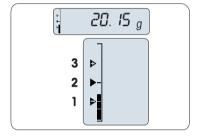

SET. I +/- %

20.28 g

#### Exit current application


To exit the current application and to return to simple weighing mode, press and hold  $\overline{\Lambda}$  (longer than 1.5 s).

# 6.4 Application statistics



The **Statistics** application allows you to generate statistics of a series of weighing values. 1 to 999 values are possible.

- 1 Call-up **APP.LIST** by pressing and holding **F**.
- 2 Select application **STAT**. by scrolling with **S**.
- 3 Activate function **STAT.** by pressing ← .

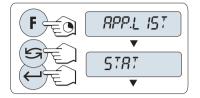

# **Memory clear question**

If the memory is already cleared (sample counter is 0) the memory clear question will not be displayed.

- 1 To continue the last statistics press ← to confirm **CLR.M:NO**.
- 2 For a new statistical evaluation clear the memory. Press to select **CLR.M:YES** and press to confirm.



#### Weighing the first sample weight


- 1 Press  $\rightarrow 0/T \leftarrow$  to zero/tare the balance if needed.
- 2 Load the first sample weight.
- 3 Press ← I. The display shows the sample count 1 and the current weight is stored as sample and the weight is printed out.

  Note: When the sample counter is displayed you may press C to undo (drop) this sample.
- 4 Unload the first sample weight.

#### Weighing further sample weights

The same procedure as for the first sample weight.

- 1...999 samples are possible.
- The next value will be accepted if the sample weight is in the range 70% – 130% of the current average value. OUT OF RANGE will be displayed if the sample is not accepted.

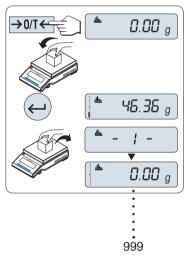

#### **Results**




- Press 🗏 if the numbers of sample are greater than or equal to 2.
- ⇒ Results are displayed and printed.

Jewelry Balances Applications

# **Displayed results**

- 1 Press ← to show the next statistical value.
- 2 Press C to cancel displaying results and to continue weighing next sample.

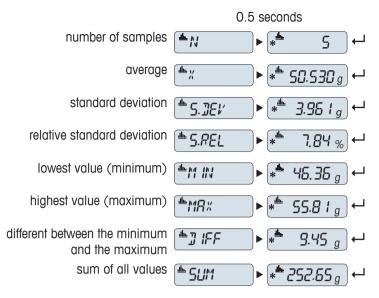

#### **Exit current application**


To exit the current application and to return to simple weighing mode, press and hold  $\Delta \bar{\Delta}$  (longer than 1.5 s).

# 6.5 Application totaling



The **TOTALING** application allows you to weigh in different samples to add their weight values and to totalize them. 1 to 999 samples are possible.


- 1 Call-up **APP.LIST** by pressing and holding **F**.
- 2 Select application **TOTAL** by scrolling with **S**.
- 3 Activate function **TOTAL** by pressing ← J.

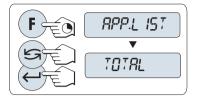

# **Memory clear question**

If the memory is already cleared (sample counter is 0) the memory clear question will not be displayed.

- 1 To continue the totaling evaluation press ← to confirm CLR.M:NO.
- 2 For a new totaling evaluation clear the memory. Press of to select **CLR.M:YES** and press to confirm.

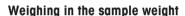

- 1 If using a container: place empty container on the weighing pan and press  $\rightarrow 0/T \leftarrow$  to zero or tare the balance.
- 2 Load the first sample weight.
- 3 Press —I. The display shows the sample count 1 and the current weight is stored.

**Note:** When the sample counter is displayed you may press **C** to undo (drop) this sample.


4 Unload the first sample weight. The display shows zero.

#### Weighing in further sample weights

The same procedure as for the first sample weight.

• 1...999 samples are possible.

# CLR.MINO CLERRE I

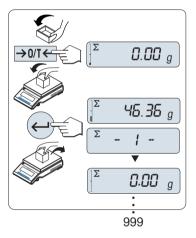

43

#### **Results**



- Press if the numbers of sample are greater than or equal to 2.
- ⇒ Results are displayed and printed.

# Displayed results

- Press ← briefly to show the totalized value.
- 2 Press C briefly to cancel.

# 0.5 seconds number of samples $\stackrel{\Sigma}{N}$ $\blacktriangleright$ $\stackrel{*}{*}$ 879 $\leftarrow$ totalized value $\stackrel{\Sigma}{TOTRL}$ $\blacktriangleright$ $\stackrel{*}{*}$ 879 $\rightarrow$ $\stackrel{\Xi}{A}$

#### **Exit current application**

To exit the current application and to return to simple weighing mode, press and hold  $\overline{\Delta}\overline{\Delta}$  (longer than 1.5 s).


Jewelry Balances Applications

# 6.6 Application multiplication factor weighing



The **Multiplication factor weighing** application allows you to multiply the weight value (in grams) by a predefined factor (result = factor \* weight) and have it calculated to a predefined number of decimal places.


- 1 Call-up **APP.LIST** by pressing and holding **F**.
- 2 Select application **FACTOR.M** by scrolling with S.
- 3 Activate function **FACTOR.M** by pressing \( \ldots \).

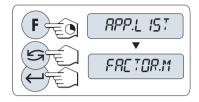

# Setting the factor value

- 1 Press to execute **SET.F.MUL**. Either the factor 1 appears as default value or the factor that was saved most recently.
- 2 To select a digit, press ← (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 3 To change the digit, press 5.
- 4 Press and hold ← to confirm the selected factor (no automatic acceptance).

**Important:** Zero for multiplication factor value is outside the allowed range, the error message **FACTOR OUT OF RANGE** will be displayed.

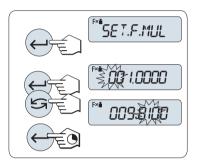


# Setting the step value

**SET.STEP** appears in the display, and the program changes automatically to allow the display increments to be entered. The smallest possible display increment appears as default value, or the last value that was saved.

- 1 Press ← to execute **SET.STEP**.
- 2 To select a digit, press ← (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 3 To change the digit, press 5.
- 4 Press and hold ← to confirm the selected step (no automatic acceptance).

**Important:** The allowed range for the step depends on the factor and the resolution of the balance. If it is outside the allowed range the error message **STEP OUT OF RANGE** will be displayed.

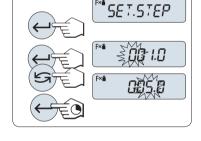

**Important:** If without any key press within 60 seconds, the balance return to the previous active application. Press **C** to cancel and returns to the previous active application.

On completion of the setting procedure, your balance is ready for multiplication factor weighing.

#### Weighing procedure


- 1 Press  $\rightarrow 0/T \leftarrow$  to zero/tare.
- 2 Load sample weight on weighing pan.
- 3 Read the result. The appropriate calculation is then made using the weight of sample and the selected factor, the result being displayed with the selected display step.

Note: No units are displayed.

4 Unload sample weight.

# Toggling between displaying the calculated value and the measured weight

You can use the key to toggle between the calculated Value, weight value **UNIT 1**, **RECALL** value (if selected) and weight value **UNIT 2** (if different from **UNIT 1**).

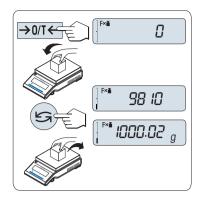

45

#### **Exit current application**

To exit the current application and to return to simple weighing mode, press and hold  $\Delta \bar{\Delta}$  (longer than 1.5 s).

Jewelry Balances Applications

# 6.7 Application division factor weighing



The **Division factor weighing** divide a predefined factor by the weight value (in grams) (result = factor / weight) and have it rounded to a predefined number of decimal places.

- Call-up **APP.LIST** by pressing and holding **F**.
- 2 Select application **FACTOR.D** by scrolling with
- 3 Activate function **FACTOR.D** by pressing ← J.



# Setting the factor value

- 1 Press to execute **SET.F.DIV**. Either the factor 1 appears as default value or the factor that was saved most recently.
- 2 To select a digit, press (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 3 To change the digit, press 5.
- 4 Press and hold ← to confirm the selected factor (no automatic acceptance).

**Important:** Zero for division factor value is outside the allowed range, the error message FACTOR OUT OF RANGE will be displayed.



SET.STEP appears in the display, and the program changes automatically to allow the display increments to be entered. The smallest possible display increment appears as default value, or the last value that was saved.


**SET.STEP** appears in the display, and the program changes automatically to allow the display increments to be entered. The smallest possible display increment appears as default value, or the last value that was saved.


- 1 Press ← to execute **SET.STEP**.
- 2 To select a digit, press ← (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 3 To change the digit, press 5.
- 4 Press ← to confirm the selected step (no automatic acceptance).


Important: The allowed range for the step depends on the factor and the resolution of the balance. If it is outside the allowed range the error message STEP OUT OF RANGE will be displayed.

**Important:** If without any key press within 60 seconds, the balance return to the previous active application. Press **C** to cancel and returns to the previous active application.

On completion of the setting procedure, your balance is ready for division factor weighing.

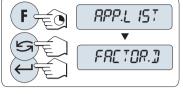

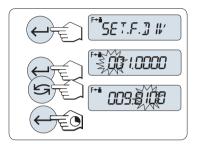

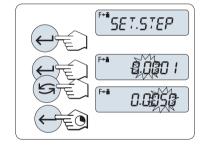

#### Weighing procedure


- 1 Press  $\rightarrow 0/T \leftarrow$  to zero/tare.
- 2 Load sample weight on weighing pan.
- 3 Read the result. The appropriate calculation is then made using the weight of sample and the selected factor, the result being displayed with the selected display step.

**Note:** No units are displayed. To avoid a division by zero, the factor division is not calculated at zero.

4 Unload sample weight.

# Toggling between displaying the calculated value and the measured weight

You can use the key to toggle between the calculated Value, weight value **UNIT 1**, **RECALL** value (if selected) and weight value **UNIT 2** (if different from **UNIT 1**).

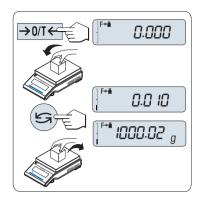

#### **Exit current application**

To exit the current application and to return to simple weighing mode, press and hold  $\vec{\Delta}$  (longer than 1.5 s).

Jewelry Balances Applications

# 6.8 Application density



The **Density** application allows you to determine the density of solid bodies and liquids. Determination of the density uses **Archimedes' principle** according to which a body immersed in a fluid undergoes an apparent loss in weight which is equal to the weight of the fluid it displaces.

To determine the density of solid bodies, we recommend you to work with the optional density kit which contains all the attachments and aids needed for convenient and precise density determination. To determine the density of liquids, you additionally need a sinker which you can also obtain from your METTLER TOLEDO dealer.


#### Note for performing of density determinations

- You can also use the hanger for weighing below the balance which belongs to your balance
- If a METTLER TOLEDO printer is attached to your balance, the settings will be automatically recorded.



We recommend you to consult the operating instructions enclosed with the density kit.

- 1 Call-up APP.LIST by pressing and holding F.
- 2 Select application **DENSITY** by scrolling with **S**.
- 3 Activate function **DENSITY** by pressing ← I.

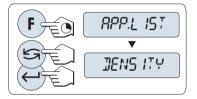

#### Setting the method for density determination

1 Select:

**SOLID**, the function for the density determination of solids, or **LIQUID**, the function for the density determination of liquids with a sinker.

2 Press ← to confirm the selection

# # SOL II

#### Exit current application


To exit the current application and to return to simple weighing mode, press and hold  $\overline{\Lambda}$  (longer than 1.5 s).

#### 6.8.1 Density determination of solids

Requirement: the method SOLID is set.

#### Setting the parameter of the auxiliary liquid

- Select the auxiliary liquid by scrolling with :: H-2-0 for distilled water, ETHANOL or FREE for a freely definable auxiliary liquid.
- 2 Press ← to confirm the selection.

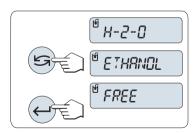

#### If you have selected water or ethanol as the auxiliary liquid

- 1 Enter the current temperature of the auxiliary liquid (read off on thermometer).
- 2 Set the value in °C. The temperature ranges from 10 °C to 30.9 °C.
- 3 To select a digit, press ← (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 4 To change the digit, press 5.
- 5 Press and hold \leftarrow to confirm the value.

**Note:** the densities of distilled water and ethanol in the range 10 °C to 30.9 °C are stored in the balance.

#### If you have selected a freely definable auxiliary liquid

- 1 Enter the density of the auxiliary liquid in g/cm<sup>3</sup> at the current temperature (read off on thermometer).
- 2 To select a digit, press  $\leftarrow$  (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 3 To change the digit, press 🥰.
- 4 Press and hold \to confirm the value.

**Important:** if without any key press within 60 seconds or by pressing **C**, the balance returns to the previous active application.

On completion of the settings, your balance is ready for performing the density determination of liquids.

Note: taring the balance is possible at any time.

The balance prompts you: PRESS ENTER TO START.

Press ← to start. Tare/Zero is executed.

The balance prompts you to weigh the solid in air WEIGH IN AIR.

- 1 Load the solid.
- 2 Press to initiate the measurement.

The balance prompts you to weigh the solid in the auxiliary liquid **WEIGH IN LIQUID**.

- 1 Load the solid.
- 2 Press ← to initiate the measurement.

The balance now shows the determined density of the solid in g/cm<sup>3</sup>.


#### Note


- This result has already been corrected for the air buoyancy. The buoyancy caused by the two immersed wires (ø 0.6 mm) can be neglected.
- By pressing C, the balance returns to PRESS ENTER TO START.

#### Result



- Press =
  - ⇒ The result will be printed.

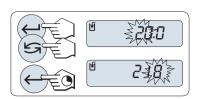

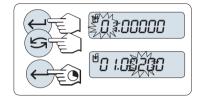









49

Jewelry Balances Applications


#### 6.8.2 Density determination of liquids

Requirement: the method LIQUID is set.

#### Setting the displacement volume of your sinker

Press and hold  $\leftarrow$  to confirm the default value of 10.0 cm<sup>3</sup> or change it if needed:

- 1 To select a digit, press ← (cyclically from left to right).
  - ⇒ The selected digit is blinking.
- 2 To change the digit, press 5.
- 3 Press and hold \to confirm the value.

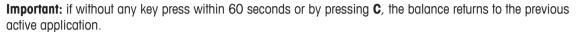

On completion of the settings, your balance is ready for performing the density determination of liquids.

Note: taring the balance is possible at any time.

The balance prompts you: PRESS ENTER TO START.

Press ← to start.



\$0.300.000

The balance prompts you to weigh the sinker in air WEIGH IN AIR.

- 1 Position the sinker.
- 2 Press ← to initiate the measurement.

The balance prompts you to weigh the sinker in the liquid **WEIGH IN LIQUID**.

- Pour the liquid into the beaker. Make sure that the sinker is immersed by al least 1 cm in the liquid, and that there are no air bubbles in the container.
- 2 Press ← to initiate the measurement.

The balance now shows the determined density of the liquid at the current temperature (read off on the thermometer).

#### Note

- This result has already been corrected for the air buoyancy. The buoyancy caused by the immersed wire (Ø 0.2 mm) of the sinker can be neglected.
- By pressing **C**, the balance returns to **PRESS ENTER TO START**.

#### Result



- Press =
  - ⇒ The result will be printed.

#### 6.8.3 Formulae used to calculate density

The **DENSITY**"application is based on the formulae listed below.









#### Formulae for determining the density of solids with compensation for air density

$$\rho = \frac{A}{A-B} (\rho_0 - \rho_L) + \rho_L$$

$$V = \alpha \frac{A - B}{\rho_0 - \rho_1}$$

 $\rho$  = Density of the sample

A = Weight of the sample in air

B = Weight of the sample in the auxiliary liquid

V = Volume of the sample

 $\rho_0 \quad = \quad \text{ Density of the auxiliary liquid}$ 

 $\rho_L$  = Density of air (0.0012 g/cm<sup>3</sup>)

 $\alpha$  = Weight correction factor (0.99985), to take the atmospheric buoyancy of the adjustment weight into account

#### Formula for determining the density of liquids with compensation for air density

$$\rho = \alpha \frac{P}{V} + \rho_L$$

 $\rho$  = Density of the liquid

Weight of the displaced liquid

V = Volume of the sinker

 $\rho_L$  = Density of air (0.0012 g/cm<sup>3</sup>)

 $\alpha$  = Weight correction factor (0.99985), to take the atmospheric buoyancy of the adjustment weight into account

T/°C	0.0	0.1	0.2	0.3	0.4	0.5	0.6	0.7	0.8	0.9
10.	0.99973	0.99972	0.99971	0.99970	0.99969	0.99968	0.99967	0.99966	0.99965	0.99964
11.	0.99963	0.99962	0.99961	0.99960	0.99959	0.99958	0.99957	0.99956	0.99955	0.99954
12.	0.99953	0.99951	0.99950	0.99949	0.99948	0.99947	0.99946	0.99944	0.99943	0.99942
13.	0.99941	0.99939	0.99938	0.99937	0.99935	0.99934	0.99933	0.99931	0.99930	0.99929
14.	0.99927	0.99926	0.99924	0.99923	0.99922	0.99920	0.99919	0.99917	0.99916	0.99914
15.	0.99913	0.99911	0.99910	0.99908	0.99907	0.99905	0.99904	0.99902	0.99900	0.99899
16.	0.99897	0.99896	0.99894	0.99892	0.99891	0.99889	0.99887	0.99885	0.99884	0.99882
17.	0.99880	0.99879	0.99877	0.99875	0.99873	0.99871	0.99870	0.99868	0.99866	0.99864
18.	0.99862	0.99860	0.99859	0.99857	0.99855	0.99853	0.99851	0.99849	0.99847	0.99845
19.	0.99843	0.99841	0.99839	0.99837	0.99835	0.99833	0.99831	0.99829	0.99827	0.99825
20.	0.99823	0.99821	0.99819	0.99817	0.99815	0.99813	0.99811	0.99808	0.99806	0.99804
21.	0.99802	0.99800	0.99798	0.99795	0.99793	0.99791	0.99789	0.99786	0.99784	0.99782
22.	0.99780	0.99777	0.99775	0.99773	0.99771	0.99768	0.99766	0.99764	0.99761	0.99759
23.	0.99756	0.99754	0.99752	0.99749	0.99747	0.99744	0.99742	0.99740	0.99737	0.99735
24	0.99732	0.99730	0.99727	0.99725	0.99722	0.99720	0.99717	0.99715	0.99712	0.99710
25.	0.99707	0.99704	0.99702	0.99699	0.99697	0.99694	0.99691	0.99689	0.99686	0.99684
26.	0.99681	0.99678	0.99676	0.99673	0.99670	0.99668	0.99665	0.99662	0.99659	0.99657
27.	0.99654	0.99651	0.99648	0.99646	0.99643	0.99640	0.99637	0.99634	0.99632	0.99629
28.	0.99626	0.99623	0.99620	0.99617	0.99614	0.99612	0.99609	0.99606	0.99603	0.99600
29.	0.99597	0.99594	0.99591	0.99588	0.99585	0.99582	0.99579	0.99576	0.99573	0.99570
30.	0.99567	0.99564	0.99561	0.99558	0,99555	0,99552	0.99549	0.99546	0.99543	0.99540

Jewelry Balances Applications

T/°C	0.0	0.1	0.2	0.3	0.4	0.5	0.6	0.7	0.8	0.9
10.	0.79784	0.79775	0.79767	0.79758	0.79750	0.79741	0.79733	0.79725	0.79716	0.79708
11.	0.79699	0.79691	0.79682	0.79674	0.79665	0.79657	0.79648	0.79640	0.79631	0.79623
12.	0.79614	0.79606	0.79598	0.79589	0.79581	0.79572	0.79564	0.79555	0.79547	0.79538
13.	0.79530	0.79521	0.79513	0.79504	0.79496	0.79487	0.79479	0.79470	0.79462	0.79453
14.	0.79445	0.79436	0.79428	0.79419	0.79411	0.79402	0.79394	0.79385	0.79377	0.79368
15.	0.79360	0.79352	0.79343	0.79335	0.79326	0.79318	0.79309	0.79301	0.79292	0.79284
16.	0.79275	0.79267	0.79258	0.79250	0.79241	0.79232	0.79224	0.79215	0.79207	0.79198
17.	0.79190	0.79181	0.79173	0.79164	0.79156	0.79147	0.79139	0.79130	0.79122	0.79113
18.	0.79105	0.79096	0.79088	0.79079	0.79071	0.79062	0.79054	0.79045	0.79037	0.79028
19.	0.79020	0.79011	0.79002	0.78994	0.78985	0.78977	0.78968	0.78960	0.78951	0.78943
20.	0.78934	0.78926	0.78917	0.78909	0.78900	0.78892	0.78883	0.78874	0.78866	0.78857
21.	0.78849	0.78840	0.78832	0.78823	0.78815	0.78806	0.78797	0.78789	0.78780	0.78772
22.	0.78763	0.78755	0.78746	0.78738	0.78729	0.78720	0.78712	0.78703	0.78695	0.78686
23.	0.78678	0.78669	0.78660	0.78652	0.78643	0.78635	0.78626	0.78618	0.78609	0.78600
24.	0.78592	0.78583	0.78575	0.78566	0.78558	0.78549	0.78540	0.78532	0.78523	0.78515
25.	0.78506	0.78497	0.78489	0.78480	0.78472	0.78463	0.78454	0.78446	0.78437	0.78429
26.	0.78420	0.78411	0.78403	0.78394	0.78386	0.78377	0.78368	0.78360	0.78351	0.78343
27.	0.78334	0.78325	0.78317	0.78308	0.78299	0.78291	0.78282	0.78274	0.78265	0.78256
28.	0.78248	0.78239	0.78230	0.78222	0.78213	0.78205	0.78196	0.78187	0.78179	0.78170
29.	0.78161	0.78153	0.78144	0.78136	0.78127	0.78118	0.78110	0.78101	0.78092	0.78084
30.	0.78075	0.78066	0.78058	0.78049	0.78040	0.78032	0.78023	0.78014	0.78006	0.77997

Density of  $\mathrm{C_2H_5OH}$  according to the "American Institute of Physics Handbook".

# 7 Communication with Peripheral Devices

#### 7.1 Function PC-Direct

The numerical value displayed at the balance can be transferred to the cursor position in Windows Applications (e.g. Excel, Word) as by typing with the keyboard.

Note: The units will not be transferred.

#### Requirements

- PC with one of the Microsoft Windows® operating system 32bit/64bit: XP (SP3), Vista (SP2), Win 7 (SP1) or Win 8.
- · Serial interface RS232 or USB.
- Administrator rights for installing software (for USB not required).
- Windows application (e.g. Excel).
- Balance to PC connection with cable RS232 or USB.

#### Settings at the balance



## **NOTICE**

Disconnect the USB connection from the balance prior to changing settings.

Balance interface settings, see Interface menu:


- Topic RS232 or USB: set PC-DIR. and select the most appropriate option for the desired weighing
  result.
- Topic RS.TX.E.O.L./RS E.O.L. or USB E.O.L./USB E.O.L:
  - set **<TAB>** to write into the same row (e.g. in Excel).
  - set <CR><LF> to write into the same column (e.g. in Excel).
- · Save changes.

#### Installing SerialPortToKeyboard

Operation of PC-Direct via serial port RS232 requires the installation of **SerialPortToKeyboard** on your host computer. The file **SerialPortToKeyboard.exe** can be found on the CD-ROM in the folder Software. If you have any questions please contact a METTLER TOLEDO representative.

#### Download SerialPortToKeyboard

- 1 Insert the product CD in the CD/DVD drive of the host computer.
- 2 Tap Software.
- 3 Download and unpack the .exe file from the CD on your computer.
- 4 Click **Save** to download to your specified location.
- 5 Right-click on the downloaded install program **SerialPort- ToKeyboard.exe** and select **Run as Administrator** from the menu.
- 6 If a safety warning appears, allow Windows to install.
- 7 Click **Next** and follow the installer's instructions.

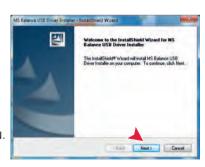


#### Settings at the PC

#### Settings for SerialPortToKeyboard

- Select the serial port COM to be used for connection with the balance.
- 2 Set the Baud Rate to 9600.
- 3 Activate Connect

#### Note

- The window can be minimized.
- Closing of the window terminates the session.

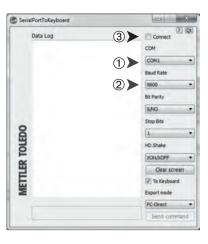

# **Checking operation**

- 1 Start SerialPortToKeyboard (RS232)
- 2 Start Excel (or another application) at the PC.
- 3 Activate a cell in Excel.

According to your selected **PC-DIR.** option, the displayed values will appear e.g. in the column one after the other one in the different rows.

# 8 Error and Status Messages

# 8.1 Error messages

Error messages in the display draw your attention to incorrect operation or that the balance could not execute a procedure properly.

Error message	Cause	Rectification
NO STABILITY	No stability.	Ensure more stable ambient conditions. If not possible, check settings for environment.
WRONG ADJUSTMENT WEIGHT	Wrong adjustment weight on pan or none at all.	Place required adjustment weight in center of pan.
REFERENCE TOO SMALL	Reference for piece counting too small.	Increase reference weight.
EEPROM ERROR - PLEASE CONTACT CUSTOMER SERVICE	<ul> <li>EEPROM (memory) error.</li> <li>Excessive mains voltage fluctuation or strong glitches occurred.</li> </ul>	Please contact METTLER TOLEDO customer service.
WRONG CELL DATA - PLEASE CONTACT CUSTOMER SERVICE	Wrong cell data.	Please contact METTLER TOLEDO customer service.
NO STANDARD ADJUSTMENT - PLEASE CONTACT CUSTOMER SERVICE	No standard calibration.	Please contact METTLER TOLEDO customer service.
PROGRAM MEMORY DEFECT - PLEASE CONTACT CUSTOMER SERVICE	Program memory defect.	Please contact METTLER TOLEDO customer service.
TEMP SENSOR DEFECT - PLEASE CONTACT CUSTOMER SERVICE	Temperature sensor defect.	Please contact METTLER TOLEDO customer service.
WRONG LOAD CELL BRAND - PLEASE CONTACT CUSTOMER SERVICE	Wrong load cell brand.	Please contact METTLER TOLEDO customer service.
WRONG TYPE DATA SET - PLEASE CONTACT CUSTOMER SERVICE	Wrong type data set.	Please contact METTLER TOLEDO customer service.
BATTERY BACKUP LOST - CHECK DATE TIME SETTINGS	Backup battery is empty. This battery ensures that the date and time are not lost when the balance is disconnected from power.	Connect the balance to the power supply for charging the battery (e.g. during the night) or contact METTLER TOLEDO customer service.
г	Overload - The weight on the pan exceeds the weighing capacity of the balance.	Reduce the weight on the weighing pan.
LJ	Underload	Check that the weighing pan is positioned correctly.
ABOVE INITIAL ZERO RANGE	Wrong weighing pan or pan is not empty.	Mount correct weighing pan or unload weighing pan.
BELOW INITIAL ZERO RANGE	Wrong weighing pan or pan is missing.	Mount correct weighing pan.
MEM.FULL	Memory full.	Clear the memory and start a new evaluation.

Jewelry Balances Error and Status Messages

Error message	Cause	Rectification
FACTOR OUT OF RANGE	Factor is outside the allow range.	Select a new factor.
STEP OUT OF RANGE	Step is outside the allowed range.	Select a new step.
	Sample weight is outside the allowed range.	Unload the pan and load a new sample weight.



# NOTICE

# Damage of the balance or software

In some countries, excessive mains voltage fluctuations and strong glitches may occur. This may affect the balance functions or damage the software.

- Use a voltage regulator for stabilizing.

# 8.2 Status messages

56

Status messages are displayed by means of small icons. The status icons indicate the following:

Status Icon	Signification
	<b>Service Reminder</b> Your balance is due for servicing. Contact your dealer's customer service department as soon as possible to have a technician service your balance. See menu topic <b>SRV.ICON</b> .

Error and Status Messages Jewelry Balances

#### 9 Maintenance

# 9.1 Cleaning and service

Every now and then, clean the weighing pan, draft shield element, bottom plate, draft shield (depending on the model) and housing of your balance. Your balance is made from high-quality, durable materials and can therefore be cleaned using a damp cloth or with a standard cleaning agent.

To thoroughly clean the draft shield glass panels, remove the draft shield from the balance. When reinstalling the draft shield, ensure that it is in the correct position.

Please observe the following notes:



#### **MARNING**

#### Danger of death or serious injury due to electric shock!

Contact with parts that contain a live current can lead to injury and death. If the balance cannot be shut down in an emergency situation, people can be injured or the balance can be damaged.

- 1 Disconnect the balance from the power supply prior to cleaning and maintenance.
- 2 Only use METTLER TOLEDO power cable, if these need to be replaced.
- 3 Make sure that no liquid enters into the balance, terminal or AC/DC adapter.
- 4 Do not open the balance, terminal or AC/DC adapter. These contain no user-serviceable parts.



# **NOTICE**

# Danger of damage to the balance due to inappropriate cleaning methods!

The balance is made from high quality, resistant materials and can be damaged by certain cleaning agents, solvents or abrasives. If liquids enter the housing they can damage the balance.

- 1 Use only water and a mild detergent to clean the balance or terminal.
- 2 Wipe off any spills immediately.
- 3 Make sure that no liquid enters the interior of the balance.

#### **Important**

Contact a METTLER TOLEDO representative to find about the service options available – regular maintenance by an authorized service engineer will ensure consistent weighing accuracy over the long term and extend the service life of the balance.

Jewelry Balances Maintenance

# 9.2 Cleaning the draft shield


It is possible to remove the sliding glass doors for cleaning or for replacing.

#### Removing or inserting sliding glass doors

- 1 Remove the handle first.
- 2 Remove the sliding glass doors.
- 3 Install the handle after insertion of the glass door.

#### **Important**

Front and rear glass panels cannot be removed.

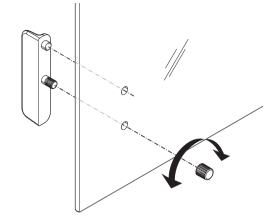

# 9.3 Disposal

58

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.



Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

Thank you for your contribution to environmental protection.

# 9.4 Firmware (Software) Updates

In the interest of its customers, METTLER TOLEDO continuously updates its internal software (firmware) for the balances. Inquire at your local METTLER TOLEDO office about upgrade and update options.

Maintenance Jewelry Balances

#### 10 Technical Data

#### 10.1 General data



# **WARNING**

### Danger of death or serious injury due to electric shock!

Contact with parts that contain a live current can lead to injury and death.

- 1 Only use an approved AC/DC adapter with a current-limited SELV output.
- 2 Ensure correct polarity ⊕ ⊕ ⊕

Standard power supply

AC adapter: Primary: 100 - 240 V, ±10%, 50/60Hz, 0.3 A

Secondary: 12 V DC, 0.84 A (with electronic overload

protection)

Balance power supply: 12 V DC, 0.84 A

Can be used up to 2000 m height above mean sea level.



# **NOTICE**

If the balance is used above 2000 m mean sea level, the optional power supply must be used.

Optional power supply

AC adapter: Primary:  $100 - 240 \text{ V, } \pm 10\%$ , 50/60 Hz

Secondary: 12 V DC ±3%, 2.5 A (with electronic overload

protection)

Cable for AC adapter: 3-core, with country-specific plug

Balance power supply: 12 V DC ±3%, 2.25 A, maximum ripple: 80 mVpp

Can be used up to 4000 m height above mean sea level.

**Protection and standards** 

Overvoltage category: II
Degree of pollution: 2

Protection: Protected against dust and water Standards for safety and EMC: See Declaration of Conformity

Range of application: For use only in closed interior rooms

**Environmental conditions** 

Height above mean sea level: Depending on the power adapter (2000 - 4000 m)

Except for China: max. 2000 m

Ambient temperature: Operating condition for ordinary lab application: +10 to 30 °C

(operability guaranteed between +5 to 40 °C)

Storage condition: -25 to 70 °C

Relative air humidity: 10% up to 80% at 31 °C, linearly decreasing to 50% at

40 °C, non-condensing

Warm-up time: At least 30 minutes (0.1 mg models 60 minutes) after

connecting the balance to the power supply

59

Jewelry Balances Technical Data

Materials

Housing: Top housing: Plastic (ABS)

Bottom housing: Die-cast aluminum, lacquered

Weighing pan: Pan ø 90 mm: Stainless steel X2CrNiMo 17-12-2 (1.4404)

All others: Stainless steel X5CrNi 18-10 (1.4301)

Draft shield element: 0.1 mg models: Stainless steel X5CrNi 18-10 (1.4301)

Draft shield: Plastic (ABS), glass

In-use-cover: Plastic (PET)

Technical Data Jewelry Balances

# 10.2 Model-specific data

# 10.2.1 Carat balances with readability of 0.001 ct / 0.1 mg

# Technical data

	JE503C	JE503CE
Limit values	1	
Maximum capacity	505 ct / 101 g	505 ct / 101 g
Readability	0.001 ct / 0.1 mg	0.001 ct / 0.1 mg
Repeatability (at nominal load)	0.1 mg	0.1 mg
Linearity deviation	0.2 mg	0.2 mg
Sensitivity temperature drift	2 ppm/°C	2 ppm/°C
Typical values		
Repeatability (at nominal load)	0.08 mg	0.08 mg
Linearity deviation	0.06 mg	0.06 mg
Minimum sample weight (U=1%, k=2)	16 mg	16 mg
Minimum sample weight OIML	10 mg	10 mg
Settling time	2 s	2 s
Adjustment	Int. Cal	Ext. Cal
Interfaces	1 RS232	1 RS232
Balance dimensions (W $\times$ D $\times$ H)	210×344×279 mm	210×344×279 mm
Weighing pan dimensions	ø 90 mm	ø 90 mm
Usable height of draft shield	170 mm	170 mm
Weight of balance	4.6 kg	4.4 kg
Weights for routine testing		
OIML CarePac	#11123002	#11123002
Weights	100 g F2, 5 g E2	100 g F2, 5 g E2
ASTM CarePac	#11123102	#11123102
Weights	100 g 1, 5 g 1	100 g 1, 5 g 1

	JE703C	JE703CE				
Limit values						
Maximum capacity	700 ct / 140 g	700 ct / 140 g				
Readability	0.001 ct / 0.1 mg	0.001 ct / 0.1 mg				
Repeatability (at nominal load)	0.1 mg	0.1 mg				
Linearity deviation	0.2 mg	0.2 mg				
Sensitivity temperature drift	2 ppm/°C	2 ppm/°C				
Typical values						
Repeatability (at nominal load)	0.08 mg	0.08 mg				
Linearity deviation	0.06 mg	0.06 mg				
Minimum sample weight (U=1%, k=2)	16 mg	16 mg				
Minimum sample weight OIML	10 mg	10 mg				
Settling time	2 s	2 s				
Adjustment	Int. Cal	Ext. Cal				
Interfaces	1 RS232	1 RS232				

61

Jewelry Balances Technical Data

	JE703C	JE703CE
Balance dimensions (W $\times$ D $\times$ H)	210×344×279 mm	210×344×279 mm
Weighing pan dimensions	ø 90 mm	ø 90 mm
Usable height of draft shield	170 mm	170 mm
Weight of balance	4.6 kg	4.4 kg
Weights for routine testing		
OIML CarePac	#11123002	#11123002
Weights	100 g F2, 5 g E2	100 g F2, 5 g E2
ASTM CarePac	#11123102	#11123102
Weights	100 g 1, 5 g 1	100 g 1, 5 g 1

	JE1103C	JE1103CE
Limit values		
Maximum capacity	1100 ct / 220 g	1100 ct / 220 g
Readability	0.001 ct / 0.1 mg	0.001 ct / 0.1 mg
Repeatability (at nominal load)	0.1 mg	0.1 mg
Linearity deviation	0.2 mg	0.2 mg
Sensitivity temperature drift	2 ppm/°C	2 ppm/°C
Typical values		
Repeatability (at nominal load)	0.08 mg	0.08 mg
Linearity deviation	0.06 mg	0.06 mg
Minimum sample weight (U=1%, k=2)	16 mg	16 mg
Minimum sample weight OIML	10 mg	10 mg
Settling time	2 s	2 s
Adjustment	Int. Cal	Ext. Cal
Interfaces	1 RS232	1 RS232
Balance dimensions (W $\times$ D $\times$ H)	210×344×279 mm	210×344×279 mm
Weighing pan dimensions	ø 90 mm	ø 90 mm
Usable height of draft shield	170 mm	170 mm
Weight of balance	4.6 kg	4.4 kg
Weights for routine testing		
OIML CarePac	#11123001	#11123001
Weights	200 g F2, 10 g F1	200 g F2, 10 g F1
ASTM CarePac	#11123001	#11123101
Weights	200 g F2, 10 g F1	200 g 1, 10 g 1

Technical Data Jewelry Balances

# 10.2.2 Gold balances with readability of 1 mg

# Technical data

	JE203G	JE203GE
Limit values		
Maximum capacity	220 g	220 g
Readability	1 mg	1 mg
Repeatability (at nominal load)	1 mg	1 mg
Linearity deviation	2 mg	2 mg
Sensitivity temperature drift	2 ppm/°C	2 ppm/°C
Typical values		
Repeatability (at nominal load)	0.7 mg	0.7 mg
Linearity deviation	0.6 mg	0.6 mg
Minimum sample weight (U=1%, k=2)	140 mg	140 mg
Minimum sample weight OIML	20 mg	20 mg
Settling time	1.5 s	1.5 s
Adjustment	Int. Cal	Ext. Cal
Interfaces	1 RS232	1 RS232
Balance dimensions (W $\times$ D $\times$ H)	210×319×289 mm	210×319×289 mm
Weighing pan dimensions	ø 120 mm	ø 120 mm
Usable height of draft shield	170 mm	170 mm
Weight of balance	4.6 kg	4.4 kg
Weights for routine testing		
OIML CarePac	#11123001	#11123001
Weights	200 g F2, 10 g F1	200 g F2, 10 g F1
ASTM CarePac	#11123101	#11123101
Weights	200 g 1, 10 g 1	200 g 1, 10 g 1

	JE303G	JE303GE
Limit values		
Maximum capacity	320 g	320 g
Readability	1 mg	1 mg
Repeatability (at nominal load)	1 mg	1 mg
Linearity deviation	2 mg	2 mg
Sensitivity temperature drift	2 ppm/°C	2 ppm/°C
Typical values		
Repeatability (at nominal load)	0.7 mg	0.7 mg
Linearity deviation	0.6 mg	0.6 mg
Minimum sample weight (U=1%, k=2)	140 mg	140 mg
Minimum sample weight OIML	20 mg	20 mg
Settling time	1.5 s	1.5 s
Adjustment	Int. Cal	Ext. Cal
Interfaces	1 RS232	1 RS232
Balance dimensions (W $\times$ D $\times$ H)	210×319×289 mm	210×319×289 mm
Weighing pan dimensions	ø 120 mm	ø 120 mm

Jewelry Balances Technical Data 63

	JE303G	JE303GE
Usable height of draft shield	170 mm	170 mm
Weight of balance	4.6 kg	4.4 kg
Weights for routine testing		
OIML CarePac	#11123001	#11123001
Weights	200 g F2, 10 g F1	200 g F2, 10 g F1
ASTM CarePac	#11123101	#11123101
Weights	200 g 1, 10 g 1	200 g 1, 10 g 1

	JE503G	JE503GE	
Limit values			
Maximum capacity	520 g	520 g	
Readability	1 mg	1 mg	
Repeatability (at nominal load)	1 mg	1 mg	
Linearity deviation	2 mg	2 mg	
Sensitivity temperature drift	3 ppm/°C	3 ppm/°C	
Typical values			
Repeatability (at nominal load)	0.7 mg	0.7 mg	
Linearity deviation	0.6 mg	0.6 mg	
Minimum sample weight (U=1%, k=2)	140 mg	140 mg	
Minimum sample weight OIML	20 mg	20 mg	
Settling time	1.5 s	1.5 s	
Adjustment	Int. Cal	Ext. Cal	
Interfaces	1 RS232	1 RS232	
Balance dimensions (W $\times$ D $\times$ H)	210×319×289 mm	210×319×289 mm	
Weighing pan dimensions	ø 120 mm	ø 120 mm	
Usable height of draft shield	170 mm	170 mm	
Weight of balance	4.6 kg	4.4 kg	
Weights for routine testing			
OIML CarePac	#11123007	#11123007	
Weights	500 g F2, 20 g F1	500 g F2, 20 g F1	
ASTM CarePac	#11123107	#11123107	
Weights	500 g 1, 20 g 1	500 g 1, 20 g 1	

Technical Data Jewelry Balances

# 10.2.3 Gold balances with readability of 10 mg

# Technical data

	JE1002G	JE1002GE	
Limit values			
Maximum capacity	1200 g	1200 g	
Readability	10 mg	10 mg	
Repeatability (at nominal load)	10 mg	10 mg	
Linearity deviation	20 mg	20 mg	
Sensitivity temperature drift	3 ppm/°C	3 ppm/°C	
Typical values			
Repeatability (at nominal load)	7 mg	7 mg	
Linearity deviation	6 mg	6 mg	
Minimum sample weight (U=1%, k=2)	1.4 g	1.4 g	
Minimum sample weight OIML	500 mg	500 mg	
Settling time	1 s	1 s	
Adjustment	Int.Cal	Ext.Cal	
Interfaces	1 RS232	1 RS232	
Balance dimensions (W $\times$ D $\times$ H)	200×319×100 mm	200×319×100 mm	
Weighing pan dimensions	180×180 mm	180×180 mm	
Weight of balance	3.8 kg	3.2 kg	
Weights for routine testing			
OIML CarePac	#11123008	#11123008	
Weights	1000 g F2, 50 g F2	1000 g F2, 50 g F2	
ASTM CarePac	#11123108	#11123108	
Weights	1000 g 1, 50 g 1	1000 g 1, 50 g 1	

	JE2002G	JE2002GE
Limit values	'	<u>'</u>
Maximum capacity	2200 g	2200 g
Readability	10 mg	10 mg
Repeatability (at nominal load)	10 mg	10 mg
Linearity deviation	20 mg	20 mg
Sensitivity temperature drift	3 ppm/°C	3 ppm/°C
Typical values		
Repeatability (at nominal load)	7 mg	7 mg
Linearity deviation	6 mg	6 mg
Minimum sample weight (U=1%, k=2)	1.4 g	1.4 g
Minimum sample weight OIML	500 mg	500 mg
Settling time	1 s	1 s
Adjustment	Int.Cal	Ext.Cal
Interfaces	1 RS232	1 RS232
Balance dimensions (W $\times$ D $\times$ H)	200×319×100 mm	200×319×100 mm
Weighing pan dimensions	180×180 mm	180×180 mm
Weight of balance	3.8 kg	3.2 kg

Jewelry Balances Technical Data

		JE2002G	JE2002GE
Weights for routine testing			
OIML CarePac		#11123009	#11123009
We	eights	2000 g F2, 100 g F2	2000 g F2, 100 g F2
ASTM CarePac		#11123109	#11123109
We	eights	2000 g 1, 100 g 1	2000 g 1, 100 g 1

	JE3002G	JE3002GE	
Limit values			
Maximum capacity	3200 g	3200 g	
Readability	10 mg	10 mg	
Repeatability (at nominal load)	10 mg	10 mg	
Linearity deviation	20 mg	20 mg	
Sensitivity temperature drift	3 ppm/°C	3 ppm/°C	
Typical values			
Repeatability (at nominal load)	7 mg	7 mg	
Linearity deviation	6 mg	6 mg	
Minimum sample weight (U=1%, k=2)	1.4 g	1.4 g	
Minimum sample weight OIML	500 mg	500 mg	
Settling time	1 s	1 s	
Adjustment	Int.Cal	Ext.Cal	
Interfaces	1 RS232	1 RS232	
Balance dimensions (W $\times$ D $\times$ H)	200×319×100 mm	200×319×100 mm	
Weighing pan dimensions	180×180 mm	180×180 mm	
Weight of balance	3.8 kg	3.2 kg	
Weights for routine testing			
OIML CarePac	#11123009	#11123009	
Weights	2000 g F2, 100 g F2	2000 g F2, 100 g F2	
ASTM CarePac	#11123109	#11123109	
Weights	2000 g 1, 100 g 1	2000 g 1, 100 g 1	

	JE4002G	JE4002GE	
Limit values			
Maximum capacity	4200 g	4200 g	
Readability	10 mg	10 mg	
Repeatability (at nominal load)	10 mg	10 mg	
Linearity deviation	20 mg	20 mg	
Sensitivity temperature drift	3 ppm/°C	3 ppm/°C	
Typical values			
Repeatability (at nominal load)	7 mg	7 mg	
Linearity deviation	6 mg	6 mg	
Minimum sample weight (U=1%, k=2)	1.4 g	1.4 g	
Minimum sample weight OIML	500 mg	500 mg	

Technical Data Jewelry Balances

	JE4002G	JE4002GE	
Settling time	1 s	1 s	
Adjustment	Int.Cal	Ext.Cal	
Interfaces	1 RS232	1 RS232	
Balance dimensions (W $\times$ D $\times$ H)	200×319×100 mm	200×319×100 mm	
Weighing pan dimensions	180×180 mm	180×180 mm	
Weight of balance	3.8 kg	3.2 kg	
Weights for routine testing			
OIML CarePac	#11123010	#11123010	
Weights	2000 g F2, 200 g F2	2000 g F2, 200 g F2	
ASTM CarePac	#11123110	#11123110	
Weights	2000 g 4, 200 g 4	2000 g 4, 200 g 4	

	JE5002G	JE5002GE	
Limit values			
Maximum capacity	5200 g	5200 g	
Readability	10 mg	10 mg	
Repeatability (at nominal load)	10 mg	10 mg	
Linearity deviation	20 mg	20 mg	
Sensitivity temperature drift	3 ppm/°C	3 ppm/°C	
Typical values			
Repeatability (at nominal load)	7 mg	7 mg	
Linearity deviation	6 mg	6 mg	
Minimum sample weight (U=1%, k=2)	1.4 g	1.4 g	
Minimum sample weight OIML	500 mg	500 mg	
Settling time	1 s	1 s	
Adjustment	Int.Cal	Ext.Cal	
Interfaces	1 RS232	1 RS232	
Balance dimensions (W $\times$ D $\times$ H)	200×319×100 mm	200×319×100 mm	
Weighing pan dimensions	180×180 mm	180×180 mm	
Weight of balance	3.8 kg	3.2 kg	
Weights for routine testing			
OIML CarePac	#11123011	#11123011	
Weights	5000 g F2, 200 g F2	5000 g F2, 200 g F2	
ASTM CarePac	#11123111	#11123111	
Weights	5000 g 4, 200 g 4	5000 g 4, 200 g 4	

Jewelry Balances Technical Data


# 10.3 Dimensions

# 10.3.1 Carat balances with readability of 0.001 ct / 0.1 mg

# Models:

JE503C JE503CE JE703C JE703CE JE1103C JE1103CE


68

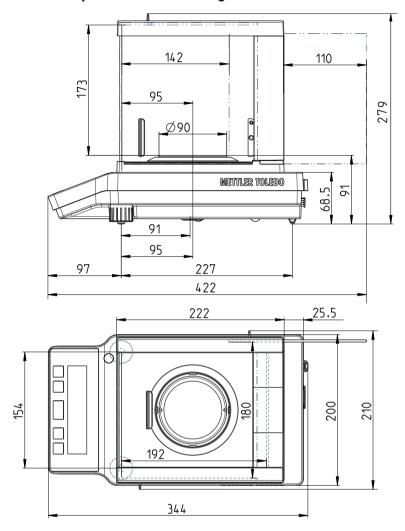

Technical Data Jewelry Balances

# 10.3.2 Gold balances with readability of 1 mg



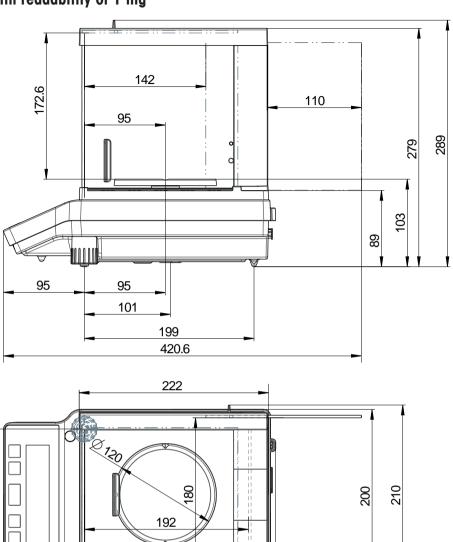



Jewelry Balances Technical Data

# 10.3.3 Gold balances with readability of 10 mg

# Models: JE1002G JE1002GE JE2002G JE2002GE JE3002G JE3002GE JE4002G JE4002G JE5002G JE5002GE

70

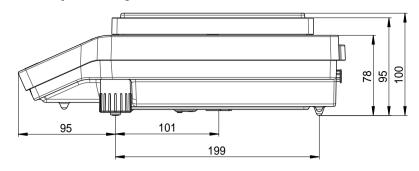

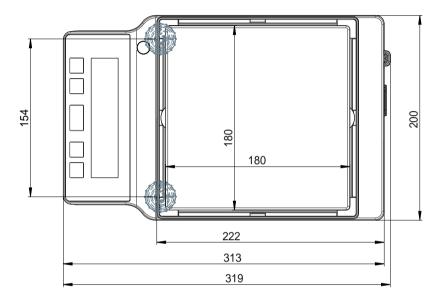

Technical Data Jewelry Balances

## 10.4 Interface specification

## 10.4.1 RS232C interface

Each balance is equipped with an RS232C Interface as standard for the attachment of a peripheral device (e.g. printer or computer).

Schematic	Item	Specification
	Interface type	Voltage interface according to EIA RS-232C/ DIN66020 CCITT V24/V.28)
DATA	Max. cable length	15 m
RXD IN TXD OUT	Signal level	Outputs: +5 V +15 V (RL = 3-7 k $\Omega$ ) -5 V15 V (RL = 3-7 k $\Omega$ ) Inputs: +3 V +25 V -3 V25 V
5 0 0 0 01	Connector	Sub-D, 9-pole, female
9 6 6 HAND SHAKE  CTS IN  RTS OUT	Operating mode	Full duplex
	Transmission mode	Bit-serial, asynchronous
	Transmission code	ASCII
	Baud rates	600, 1200, 2400, 4800, 9600, 19200, 38400 (software selectable)
POWER	Bits/parity	7-bit/none, 7-bit/even, 7-bit/odd, 8-bit/none (software selectable)
+12V OUT	Stop bits	1 stop bit
2nd display mode only	Handshake	None, XON/XOFF, RTS/CTS (software selectable)
	End-of-line	<cr><lf>, <cr>, <lf> (software selectable)</lf></cr></lf></cr>
	Power supply for 2nd display	+ 12 V, max 40 mA (software selectable, 2nd display mode only)

## 10.4.2 MT-SICS interface commands and functions

Many of the instruments and balances used have to be capable of integration in a complex computer or data acquisition system.

To enable you to integrate balances in your system in a simple manner and utilize their capabilities to the full, most balance functions are also available as appropriate commands via the data interface.

All new METTLER TOLEDO balances launched on the market support the standardized command set "METTLER TOLEDO Standard Interface Command Set" (MT-SICS). The commands available depend on the functionality of the balance.

For further information please contact your METTLER TOLEDO representative.

For further information please refer to the Reference Manual MT-SICS downloadable from the Internet under

www.mt.com/jewelry

Jewelry Balances Technical Data

# 11 Accessories and Spare Parts

## 11.1 Accessories

## **Accessories**


## **Printers**

	RS-P25 printer with RS232 connection to instrument Paper roll (length: 20 m), set of 5 pcs Paper roll (length: 13 m), self-adhesive, set of 3 pcs Ribbon cartridge, black, set of 2 pcs	11124300 00072456 11600388 00065975
	RS-P26 printer with RS232 connection to instrument (with date and time)  Paper roll (length: 20 m), set of 5 pcs  Paper roll, self-adhesive (length: 13 m), set of 3 pcs	11124303 00072456 11600388
	Ribbon cartridge, black, set of 2 pcs	00065975
	RS-P28 printer with RS232 connection to instrument (with date, time and applications)	11124304
	Paper roll (length: 20 m), set of 5 pcs Paper roll, self-adhesive (length: 13 m), set of 3 pcs	00072456 11600388
	Ribbon cartridge, black, set of 2 pcs	00065975
	P-52RUE dot matrix printer with USB, RS232, and ethernet connections	30237920
	Paper roll, set of 5 pcs Paper roll, self-adhesive, set of 3 pcs Ribbon cartridge, black, set of 2 pcs	00072456 11600388 00065975
	P-56RUE thermal printer with RS232, USB and ethernet connections, simple printouts, date and time, label printing (limited).	30094673
	Paper roll, white (length: 27 m), set of 10 pcs Paper roll, white, self-adhesive (length: 13 m), set of 10 pcs	30094723 30094724
	Paper roll, white, self-adhesive labels (550 labels), set of 6 pcs	30094725
	P-58RUE thermal printer with RS232, USB and ethernet connections, simple printouts, date and time, label printing, balance applications: statistics, formulation, totaling,	30094674
6	Paper roll, white (length: 27 m), set of 10 pcs Paper roll, white, self-adhesive (length: 13 m), set of 10 pcs	30094723 30094724
	Paper roll, white, self-adhesive labels (550 labels), set of 6 pcs	30094725

Accessories and Spare Parts

Jewelry Balances

## Cables for RS232 interface

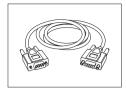


RS9 - RS9 (m/f): connection cable for PC, length = 1 m

11101051



RS9 - RS25 (m/f): connection cable for PC, length = 2 m

11101052

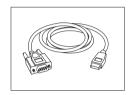

RS232 - USB converter cable — Cable with converter to connect a balance (RS232) to a USB port


64088427

## **Density determination**




Density kit ME-DNY-4 for balances with readability of 0.01 mg / 0.1 mg

30029886

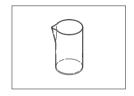

Glass beaker, height 100 mm, ø 60 mm

00238167

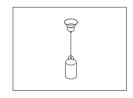

Sinker for density of liquids in conjunction with density kit Calibrated (sinker + certificate)

Recalibrated (new certificate)

00210260 00210672 00210674



Calibrated thermometer with certificate

11132685

## **Carat pans**




Carat pan XS, ø 50 mm / heigh 20 mm (set of 10 units)

12102565




Carat pan S, ø 80 mm / heigh 20 mm (set of 10 units)

12102645

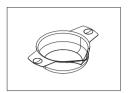

Carat pan M, ø 90 mm / heigh 30 mm (set of 10 units)

12102646

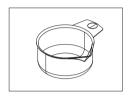

Carat pan L, ø 90 mm / heigh 45 mm (set of 10 units)

12102647

## Weighing pans



Set of weighing pan ø 160 mm with pan support for balances with readability of 10 mg and 100 mg using draft shield

30042896

#### **Draft shields**




Draft shield low with sliding doors, usable heigh 170 mm.

30042884

- for balances 0.1 mg or 1 mg
- for balances 10 mg or 100 mg, weighing pan ø 160 mm is needed (#30046407)


## **Auxiliary displays**

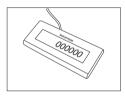

RS232 auxiliary display AD-RS-J7


12122380

#### **Protective covers**

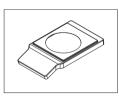

Protective cover for models with readability of 0.01 mg / 0.1 mg

30037742

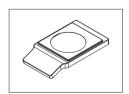


Protective cover for models with readability of 1 mg ... 100 mg

30042890


Accessories and Spare Parts Jewelry Balances

#### **Dust covers**

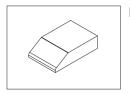

Dust cover for models without draft shield

30029051

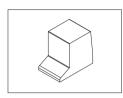

Dust cover for models with draft shield low (170 mm)

30029050


## **Anti-theft devices**



Anti-theft steel cable

11600361

## Software

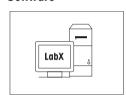

LabX direct balance (simple data transfer)

11120340

## **Transport cases**




Transport case for models with draft shield (170 mm) Transport case for models without draft shield 30046405 30046406

## Miscellaneous




Tool kit, contains brush, tweezer and glove

30046403

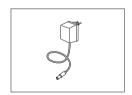


AC/DC universal adapter (EU, USA, AU, UK) 100–240 VAC, 50/60 Hz, 0.5 A, 12 VDC 1 A

11120270

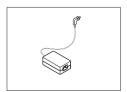

AC/DC adapter (without power cable) 100–240 V AC, 0.8 A, 50/60 Hz, 12 V DC 2.5 A

11107909

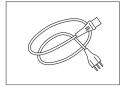

Country-specific 3-Pin power cable with grounding conductor.

Power cable AU	00088751
Power cable BR	30015268
Power cable CH	00087920
Power cable CN	30047293
Power cable DK	00087452
Power cable EU	00087925
Power cable GB	00089405
Power cable IL	00225297
Power cable IN	11600569
Power cable IT	00087457
Power cable JP	11107881
Power cable TH, PE	11107880
Power cable US	00088668
Power cable ZA	00089728

## **Adjustment weights**




OIML / ASTM Weights (with calibration certificate) see http://www.mt.com/weights

## 11.2 Spare parts

76

## JE balances with readability of 0.1 mg / 1 mg

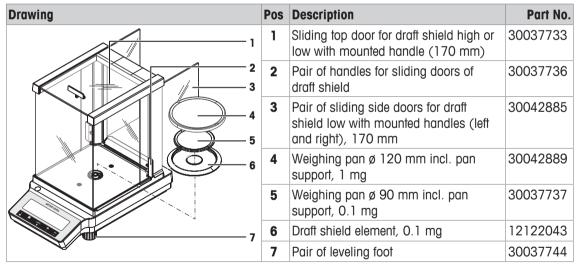

Accessories and Spare Parts

Jewelry Balances

## JE balances with readability of 10 mg

Drawing	Pos	Description	Part No.
	1	Weighing pan $180 \times 180$ mm with pan support, 10 mg	30042895
1	2	Draft shield element $180 \times 180$ mm, $10$ mg	30042897
2	3	Pair of leveling foot	30037744

## Miscellaneous

Drawing	Pos	Description	Part No.
		Bottom plate for draft shield	30037739
		Bottom plate for models without draft shield	30042901

Accessories and Spare Parts

Jewelry Balances

78

## Index

A         Data communication format         3 communication         4 communication         2 communication         4 communication         2 communication         4 communication         4 communication         2 communication         4 communication         2 communication         4 communication         3 communication         4 communication         3 communication         4 communication	Numerics		D	
Codopter	1/10 d	26	Data communication format	30
AC adapter         59         Date format         22           Accessories         72         Delivery inspection         12           Adjusting         19         Density         44           Adjustment         21         Density kit         48           Adjustment lock         28         Dimensions         68           Advanced menu         25         Disploy         39           Application check weighing         39         Disploy ponel         26           Application density         48         Dispoor ponel         26           Application icons         9         Division factor weighing         56           Application multiplication factor weighing         44         Dispooral         55           Application percent weighing         44         Proft shield         57         56           Application piece counting         41         End of line         32         40         32         40	A		Date	17, 26
Accessories         72         Delivery inspection         12           Adjusting         19         Density kit         48           Adjustment         21         Density kit         48           Adjustment lock         28         Density table for ethanol         55           Adjustment lock         28         Disploy         30           Application develweighing         39         Disploy ponel         56           Application density         48         Disploy ponel         56           Application icons         9         Disploy ponel         56           Application multiplication factor weighing         44         Draft shield         57, 56           Application percent weighing         37         Port of shield         57, 56           Application percent weighing         44         End of line         32           Application problec counting         41         End of line         32           Application statistics         41         End of line         32           Application totaling         43         Environmental conditions         55           Auto print         30         Error messages         56           Auto print         30         Error messages         56<		50	Date format	28
Adjusting 19 Density kit 44 Adjustment 21 Density kit 41 Adjustment lock 28 Dimensions 66 Advanced menu 25 Display panel 25 Application division factor weighing 39 Display panel 35 Application division factor weighing 46 Display panel 35 Application division factor weighing 46 Application division factor weighing 47 Application division factor weighing 47 Application sens 49 Division factor weighing 47 Application piece counting 48 Application piece counting 49 Application piece counting 49 Application statistics 41 Application balling 43 Application totaling 43 Application totaling 43 Application statistics 41 Application balling 43 Application balling 43 Application balling 43 Application statistics 41 Application balling 43 Application statistics 41 Application balling 43 Application balling 44 Application balling 44 Application piece counting 45 Applicati	·		Delivery inspection	12
Adjustment         21         Density tible for ethanol         5.4           Adjustment lock         28         Density toble for ethanol         5.6           Adjustment lock         28         Dimensions         68           Advanced menu         25         Disploy         33           Application check weighing         39         Disploy ponel         35           Application division factor weighing         46         Disploy ponel         56           Application in weighing         44         Disploy ponel         56           Application multiplication factor weighing         44         Draft shield         57, 56           Application percent weighing         37         Part shield         57, 56           Application statistics         41         End of line         32           Autoratic standby         28         Ethanol         52           Autoratic standby         28         Ethanol			Density	48
Adjustment lock         28         Density locite for effendol         56           Advanced menu         25         Disploy         36           Application check weighing         39         Disploy ponel         26           Application density         48         Disploy ponel         56           Application division factor weighing         46         Disposal         58           Application in untiliplication factor weighing         44         Disposal         58           Application percent weighing         37         Application percent weighing         47         57         58           Application percent weighing         34         E         6         57         58         58         58         58         59         58         59			Density kit	48
Advanced menu         25 billplay         Display         38 billplay           Application check weighing         39 bisplay increment         26 bisplay panel         27 bisplay panel         27 bisplay panel         28 bisplay panel         29 bisplay panel         28 bisplay panel         29 bis	-		Density table for ethanol	52
Application check weighing 39 Display increment 25c Application density 48 Display increment 25c Application density 48 Display ponel 25c Application in white places of the properties of the p	-		Dimensions	68
Application density			Display	30
Application division factor weighing         46         Display parily         58           Application icons         9         Division factor weighing         44           Application percent weighing         37         Port shield         57, 58           Application piece counting         34         E           Application statistics         41         End of line         32           Application statistics         41         End of line         32           Application totaling         43         Environment         27           Applications         10         Environmental conditions         55           Auto print         30         Error messages         55           Automatic standby         28         Ethanol         52           Automatic zero setting         29         External weight         20           Autozero         29         F         Punction PC-Direct         55           Basic menu         25         General safety information         44           Basic principles for operation         10         Handles         56           Bil/Parily         31         Handles         56           Bil/Parily         31         Handles         36			Display increment	26
Application icons         9 Disposition         Disposition of Section and Papel Carlos and Papel Car			Display panel	9
Application multiplication factor weighing Application percent weighing Application percent weighing Application piece counting Application piece counting Application piece counting Application statistics Application statistics Application totaling Application piece counting Application statistics  10 Environmental conditions 52 Autoprint 30 Error messages 55 Automatic standby Automatic zero setting 29 External weight 20 Autorero 29 External weight 20 Average (Statistics)  Backlight Backlight Basic menu 25 Backlight Basic menu 25 Backlight Basic menu 25 Backlight Below-the-balance weighing Bit/Parity 31 Handles Bit/Parity 31 Handles  6 Cancel 11 Char set Cancel 11 Char set Char set Cleaning 11 Char set Cleaning 11 Char set Cleaning 11 Char set Cleaning 11 Components 12 Components 12 Components 12			Disposal	58
Application multiplication factor weighing         44         Draft shield         57, 56           Application piece counting         37         E           Application statistics         41         End of line         33           Application statistics         41         End of line         33           Applications         10         Environmental conditions         55           Auto print         30         Error messages         56           Automatic standby         28         Ethanol         52           Automatic zero setting         29         External weight         20           Autoarea         29         External weight         20           Average (Statistics)         41         Factor weighing         44, 46           Backlight         28         Backlight Portor         53           Backlight principles for operation         10         Beneral safety information         4           Baudrate         31         H         H           Below-the-balance weighing         23         Handles         56           Bit/Parity         31         Handles         56           Bit/Parity         31         Handles         30           Cancel         11 <td>• •</td> <td></td> <td>·</td> <td>46</td>	• •		·	46
Application percent weighing         37           Application precent weighing         34           Application statistics         41         End of line         32           Application totaling         43         Environment         27           Applications         10         Environmental conditions         58           Auto print         30         Error messages         55           Automatic standby         28         Ethanol         52           Automatic zero setting         29         External weight         20           Autozero         29         F         Cotro weighing         44           Average (Statistics)         41         Factor weighing         44, 46           Basic menu         25         General safety information         48           Basic principles for operation         10         General safety information         48           Below-the-balance weighing         23         Handles         58           Bit/Parity         31         Handles         58           Bit/Parity         31         Handles         32           Cancel         11         Intercept the deeder         30           Change settings         11         Intercept the deeder				57, 58
Application statistics			E	
Application totaling         43         Environment         27           Applications         10         Environmental conditions         58           Auto print         30         Error messages         58           Automatic standby         28         Ethanol         52           Automatic zero setting         29         External weight         20           Autozero         29         F         Factor weighing         44, 46           B         Function PC-Direct         53           Backlight         28         General safety information         4           Basic menu         25         General safety information         4           Baudrate         31         H           Below-the-balance weighing         23         Handles         58           Bit/Parity         31         Handles         32           Concel         11         Host         32           Change settings         11         I           Char set         32         Icons         32           Check weighing         39         Identification         29           Cleaning         57         In use cover         15           Closing the menu         11		34		
Applications         10         Environmental conditions         55           Auto print         30         Error messages         55           Automatic standby         28         Ethanol         52           Automatic zero setting         29         External weight         20           Autozero         29         Factor weighing         44, 46           B         Foactor weighing         44, 46           Backlight         28         General safety information         53           Basic menu         25         General safety information         44           Below-the-balance weighing         23         Handles         58           Bit/Parity         31         Handshake         32           C         Header         30           Cancel         11         Host         30           Change settings         11         I           Char set         32         Icons         58           Check weighing         39         Identification         27           Cleaning         57         In use cover         15           Cloosing the menu         11         see protective cover         15           Components         8 <t< td=""><td>• •</td><td></td><td>End of line</td><td></td></t<>	• •		End of line	
Auto print       30       Error messages       55         Automatic standby       28       Ethanol       52         Automatic zero setting       29       External weight       20         Autozero       29       Factor weighing       44, 46         B       Founction PC-Direct       53         Backlight       28       General safety information       44         Basic menu       25       General safety information       4         Baudrate       31       H       4         Below-the-balance weighing       23       Handles       56         Bit/Parity       31       Handshake       32         C       Header       30         Cancel       11       Host       30         Change settings       11       I         Char set       32       Icons       56         Check weighing       39       Identification       27         Cleaning       57       In use cover       15         Closing the menu       11       see protective cover       15         Components       8       Input principle       11         Components       12       Components       13	Application totaling	43	Environment	
Automatic standby       28       Elhanol       52         Automatic zero setting       29       External weight       20         Autozero       29       Factor weighing       44, 46         B       Function PC-Direct       53         Backlight       28       General safety information       44         Basic menu       25       General safety information       4         Baudrate       31       H       H         Below-the-balance weighing       23       Handles       58         Bit/Parity       31       Handshake       32         C       Header       30         Cancel       11       I         Change settings       11       I         Char set       32       Icons       30         Check weighing       39       Identification       27         Cleaning       57       In use cover       15         Closing the menu       11       see protective cover       18         Components       8       Input principle       11         Conventions and symbols       3       Installing         Customer fine adjustment       21       Components       13 <td>Applications</td> <td>10</td> <td>Environmental conditions</td> <td>59</td>	Applications	10	Environmental conditions	59
Automatic zero setting       29       External weight       20         Autozero       29       F         Average (Statistics)       41       Factor weighing       44, 46         B       Function PC-Direct       53         Backlight       28       G         Basic menu       25       General safety information       4         Basic principles for operation       10       H         Budrate       31       H         Below-the-balance weighing       23       Handles       58         Bit/Parity       31       Handshake       32         Cancel       11       Host       30         Change settings       11       I         Change settings       11       I         Check weighing       39       Identification       27         Cleaning       57       In use cover         Closing the menu       11       see protective cover       18         Components       8       Input principle       11         Conventions and symbols       3       Installing         Customer fine adjustment       21       Components       13	Auto print	30	Error messages	55
Autozero         29 Average (Statistics)         #         #         Feator weighing         44, 46, 46         #         #         #         Feator weighing         44, 46         #         #         #         Feator weighing         44, 46         #         #         #         Feator weighing         Feator weighing         #         Feator weighing and safety information         #	Automatic standby	28	Ethanol	52
Average (Statistics)         41         Factor weighing         44.4 46           B         Function PC-Direct         53           Backlight         28         G           Basic menu         25         General safety information         4           Basic principles for operation         10         H           Baudrate         31         H           Below-the-balance weighing         23         Handles         58           Bit/Parity         31         Handshake         32           Cancel         11         L         L           Change settings         11         I         L           Char set         32         Icons         36           Check weighing         39         Identification         27           Cleaning         57         In use cover           Closing the menu         11         see protective cover         18           Components         8         Input principle         11           Conventions and symbols         3         Installing           Customer fine adjustment         21         Components         13	Automatic zero setting	29	External weight	20
Best (a)         Factor weighing         44, 46           Backlight         28         General safety information         4           Basic menu         25         General safety information         4           Basic principles for operation         10         H           Baudrate         31         H           Below-the-balance weighing         23         Handles         58           Bit/Parity         31         Handshake         32           Cancel         11         Host         30           Change settings         11         I           Chars set         32         Icons         9           Check weighing         39         Identification         27           Cleaning         57         In use cover         15           Closing the menu         11         see protective cover         15           Components         8         Input principle         11           Conventions and symbols         3         Installing           Customer fine adjustment         21         Components         13	Autozero	29	F	
B         Function PC-Direct         53           Backlight         28         6           Basic menu         25         General safety information         4           Basic principles for operation         10         H           Baudrate         31         H           Below-the-balance weighing         23         Handles         58           Bit/Parity         31         Handshake         32           Concel         11         Host         30           Change settings         11         I           Char set         32         Icons         9           Check weighing         39         Identification         27           Cleaning         57         In use cover         15           Closing the menu         11         see protective cover         15           Components         8         Input principle         11           Conventions and symbols         3         Installing           Customer fine adjustment         21         Components         13	Average (Statistics)	41		11 16
Backlight         28         G           Basic menu         25         General safety information         4           Basic principles for operation         10         H           Baudrate         31         H           Below-the-balance weighing         23         Handles         58           Bit/Parity         31         Handshake         32           Cancel         Header         30           Cancel         11         I           Change settings         11         I           Char set         32         Icons         9           Check weighing         39         Identification         27           Cleaning         57         In use cover         15           Closing the menu         11         see protective cover         15           Conventions and symbols         3         Installing           Customer fine adjustment         21         Components         13	В			
Basic menu         25         General safety information         4           Basic principles for operation         10         H           Baudrate         31         H           Below-the-balance weighing         23         Handles         58           Bit/Parity         31         Handshake         32           Cancel         11         Host         30           Change settings         11         I           Char set         32         Icons         9           Check weighing         39         Identification         27           Cleaning         57         In use cover         15           Components         8         Input principle         11           Conventions and symbols         3         Installing         12           Customer fine adjustment         21         Components         13	Backlight	28		00
Basic principles for operation Baudrate Below-the-balance weighing Bit/Parity Below-the-balance weighing Below-	•		<b>G</b>	
Baudrate31HBelow-the-balance weighing23Handles58Bit/Parity31Handshake32CancelHeader30Change settings11IChar set32Icons9Check weighing39Identification27Cleaning57In use coverClosing the menu11see protective cover18Components8Input principle11Conventions and symbols3InstallingCustomer fine adjustment21Components13			General safety information	4
Bit/Parity 31 Handshake 32  Cancel Header 30 Change settings 11 I Char set 32 Icons Check weighing 39 Identification 27 Cleaning 57 In use cover Closing the menu 11 see protective cover 15 Components 8 Input principle 11 Conventions and symbols 3 Installing Customer fine adjustment 21 Components 13		31	Н	
Cancel Header Host 30 Change settings 11 I Char set 32 Icons Check weighing 39 Identification 27 Cleaning 57 In use cover Closing the menu 11 see protective cover 15 Conventions and symbols 3 Installing Customer fine adjustment 21 Components 13 Components 14 Components 15 Component	Below-the-balance weighing	23	Handles	58
Cancel 11 Host 30 Change settings 11 L Char set 32 Icons 9 Check weighing 39 Identification 27 Cleaning 57 In use cover Closing the menu 11 see protective cover 15 Components 8 Input principle 11 Conventions and symbols 3 Installing Customer fine adjustment 21 Components 13	Bit/Parity	31	Handshake	32
Change settings  Char set  Char set  Check weighing  Cleaning  Cleaning  Closing the menu  Components  Conventions and symbols  Customer fine adjustment  Char set  32	C		Header	30
Change settings11Image: Change settingsChar set32IconsCheck weighing39IdentificationCleaning57In use coverClosing the menu11see protective coverComponents8Input principleConventions and symbols3InstallingCustomer fine adjustment21Components	Cancel	11	Host	30
Char set32Icons32Check weighing39Identification27Cleaning57In use coverClosing the menu11see protective cover15Components8Input principle11Conventions and symbols3InstallingCustomer fine adjustment21Components13			1	
Check weighing39Identification27Cleaning57In use coverClosing the menu11see protective cover15Components8Input principle11Conventions and symbols3InstallingCustomer fine adjustment21Components13	•	32	lcons	ç
Cleaning 57 In use cover  Closing the menu 11 see protective cover 15  Components 8 Input principle 11  Conventions and symbols 3 Installing  Customer fine adjustment 21 Components 13	Check weighing	39		
Closing the menu 11 see protective cover 15 Components 8 Input principle 11 Conventions and symbols 3 Installing Customer fine adjustment 21 Components 13	Cleaning	57		
Components8Input principle11Conventions and symbols3InstallingCustomer fine adjustment21Components13	Closing the menu	11		1.5
Conventions and symbols 3 Installing Customer fine adjustment 21 Components 13	Components	8	•	
Customer fine adjustment 21 Components 13	Conventions and symbols	3		
5 Components	Customer fine adjustment	21	•	1.3
	-		·	

Jewelry Balances Index 79

Intended use	4	Power cable	16
Interface		Power supply	59
MT-SICS	71	Power supply voltage	16
Interface menu	25, 29	Print	23
Interface RS232C	29, 71	Print menu	27
Internal weight	19	Printer	29
Interval	32	Protect	26
Introduction	3	Protection and standards	59
L		Protective cover	15
Level indicator	17	R	
Leveling	17	Recall	23, 28
Line feed	30	Reset	27
Liquid	48	RS232C interface	29, 71
Liquids	50	S	
Local conditions	16		
Location	16	Safety Information	4
Lock adjustment	28	General Intended use	4
М			4
	26	Signal words	
Manual adjustment with external weight	20	Warning Symbols	4 11
Manual adjustment with external weight Manual adjustment with internal weight	19	Saving settings Select menu	10
Materials			10
Menu	60 26	Select menu topic	10
Menu Advanced	25 25	Selecting weighing application Service	29, 57
Menu Basic	25 25	Service date reset	29, 37
Menu Interface	25, 29	Service dule reser	29
	25, 29	Service reminder	29
Menu operation  Menu protection	26	Set identification	2 <del>9</del> 27
Menu topic	11, 26	Setting Date and Time	17
MT-SICS	71	Setting bale and Time  Setting up the balance	17
Multiplication factor weighing	44	Signal words	4
	44	Signature line	30
N		Sinker	48, 50
Numerical values	11	Solids	48
0		Spare parts	77
ON / OFF	17, 23	Standard deviation (Statistics)	41
Operating temperature	17	Standby	28
Overview	8	Statistics	41
P		Status icons	9
	20	Status messages	56
PC-DIR	29 53	Stop bit	32
PC-Direct	53 27	Submenu	11
Percent weighing	37	Switching	
Performing a simple weighing	22	On	17
Piece counting	34	Switching off	23

80 Index Jewelry Balances

Switching weight units	22
Symbols	3
Symbols and conventions	3
T	
Taring	22
Technical data	59
Technical data model-specific	61
Time	17, 26
Time format	28
Topic	11, 26
Totaling	43
Transmit data	23
Transport over long distances	24
Transport over short distances	24
Transporting the balance	24
U	
Unit	26, 27
Unpacking	12
W	
Warm-up time	17
Warning Symbols	4
Weighing application	10
Weighing-in aid	23
Weight unit	22, 26, 27
Z	
Zero print	30
Zero setting	22
Zeroing	29

Jewelry Balances Index 81

82 Index Jewelry Balances



Good Weighing Practice<sup>™</sup>

GWP® is the global weighing standard, ensuring consistent accuracy of weighing processes, applicable to all equipment from any manufacturer It helps to:

- Choose the appropriate balance or scale
- Calibrate and operate your weighing equipment with security
- Comply with quality and compliance standards in laboratory and manufacturing

www.mt.com/GWP

www.mt.com/jewelry	www.m	it.com/	iewe	lrv
--------------------	-------	---------	------	-----

For more information

Mettler-Toledo GmbH

Im Langacher 44 8606 Greifensee, Switzerland www.mt.com/contact

Subject to technical changes.

© Mettler-Toledo GmbH 02/2017 30066320C en



30066320