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Welcome to the Found New Hardware Wizard		
This wizard helps you install software for: USB CDC serial port emulation		
If your hardware came with an installation CD or floppy disk, insert it now.		
What do you want the wizard to do?		
 Install from a list or specific location (Advanced) 		
Click Next to continue.		
	Welcome to the Found New Hardware Wizard This wizard helps you install software for: USB CDC serial port emulation If your hardware came with an installation CD or floppy disk, insert it now. What do you want the wizard to do? Install the software automatically (Recommended) Install from a list or specific location (Advanced) Click Next to continue.	Welcome to the Found New Hardware Wizard This wizard helps you install software for: USB CDC serial port emulation Image: Serial port emulation

Har dwa	re Installation	
<u>.</u>	The software you are installing for this hardware: USB CDC serial port emulation has not passed Windows Logo testing to verify its compatibility with Windows XP. (<u>Tell me why this testing is important.</u>) Continuing your installation of this software may impair or destabilize the correct operation of your system either immediately or in the future. Microsoft strongly recommends that you stop this installation now and contact the hardware vendor for software that has passed Windows Logo testing. Continue Anyway STOP Installation	第二步: 点击继续:

Found New Hardware Wizard	
Completing the Found New Hardware Wizard The wizard has finished installing the software for: USB CDC serial port emulation Click Finish to close the wizard.	第三步: 点击安装完毕

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