
# **Operating Instructions**

# **LC-PVolume**

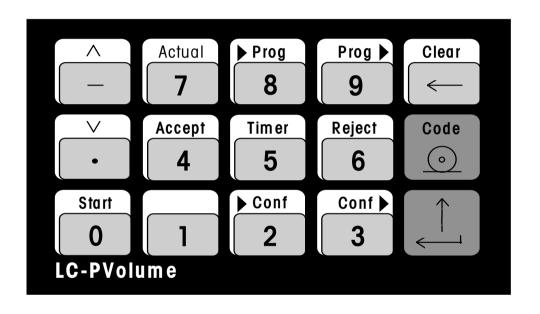

Con	tents	Page
1	METTLER TOLEDO LC-PVolume	3
2	First steps	4
	Setting up the system	4
	Checking the system using a test weight	4
3	General key functions/entries	5
4	Configuring the system	6
5	Defining the program (test parameter)	7
	Printing out program catalog	9
	Printing out test parameters	9
	Deleting test parameters	9
	Clearing individual string parameters	9
_	Default position: -59-	9
6	How does LC-PVolume work?	10
_	The Gravimetric Analysis	11
7	Error messages during operation of LC-PVolume	12
8	What if?	12
9	Sampling operation (examples)	13
10	System startup and maintenance	20
Appe	ndix A: Technical data of LC-PVolume Controller	22
Appe	ndix B: Standard and optional equipment	22
Index	(	23

#### **METTLER TOLEDO LC-PVolume**

LC-PVolume is an aid for evaluating performance of volumetric equipment. It fulfills the requirements of

- the primary test method of Guideline I8 ("Determining the Performance of Volumetric Equipment"), proposed by the National Committee for Clinical Laboratory Standards (NCCLS).
- ISO / DIS 8655/1
- DIN 12650/6

The test method, called the Gravimetric Analysis (→ section 6), is a weighing procedure. Aside from its repetitive nature, it requires a number of physical and statistical calculations, making the entire procedure a tedious one. With LC-PVolume, the length of time it takes to finish a complete test cycle would be greatly reduced. This is mainly because the operator is relieved of several tasks, namely: writing down each weighing result, calculation of results, and creating a record of the test procedure. Moreover, LC-PVolume allows for internal storage of definitions for test parameters, which further shortens the test cycle time.

Special features of the system:

Print mode

Max. number of programs 60; storage positions 0...59 (with -59- as

user specified default for defining other

empty positions)

Input: g or mg...Output: mg ( $\rightarrow \mu l$ ) Units Sample size

4/10/30

Two level password protection System configuration and programs

> Line printer, selectable in the configuration menu or via special keypress (press and hold down [1] for a few seconds while in standby mode to go to printer mode; to return to LC-PVolume mode, press and hold

down [Actual] for a few seconds).


Any one of the following METTLER TOLEDO balances may be attached:

- AM balances with standard software V10.45.XX or higher
- AT, MT, UMT balances with software version 1.09 or higher

Also for balances with restricted display function: (DeltaTrac for position and test indicator not present)

- AJ balances with data interface Option 018
- AE (Option 012) balances
- B and G series (CAN/SICS balances)
- B-S and G-S series (RS232)

#### LC-PVolume keypad

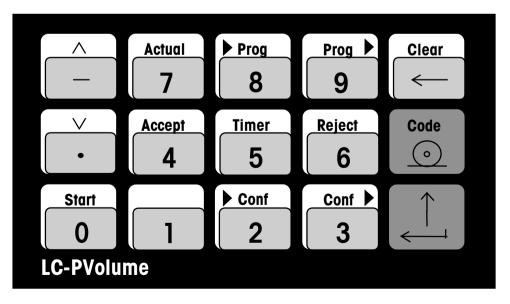

## 2 First steps

#### Setting up the system

	Unpack and position the LC-PVolume Controller					
	Install batteries ( $\rightarrow$ section 10: System startup and maintenance)					
	Insert paper (→ section 10: System startup and maintenance)					
	Connect LC-PVolume Controller and balance with the required cable					
	and attach to power supply ( $\rightarrow$ Appendix B)					
	Switch on balance and set to default settings: RESET. For some					
	certified M balances/scales, Au (Auxiliary) must also be set to ON in					
	the I-FACE menu to enable all functions (→ balance operating inst.).					
	Switch on LC-PVolume Controller and balance:					
	Status display of the LC-PVolume Controller (LED):					
	LED flashes rapidly Initialization after switching on					
	LED flashes slowly Balance is switched off, or LC-PVolume is in					
	printer mode, or there is no connection to					
	balance/scale					
	LED constantly lit up RS232C interface in order and LC-PVolume					
	, .					
П	system ready for operation  Configuration ( continued to configuring the eventum)					
ш	Configuration (→ section 4: Configuring the system)					
	Important: Enter date and time					
_	Finally, check the system using the routine opposite.					
	Define default program position -59- first ( $\rightarrow$ section 5) to have your					
	company's specific data preset on all other positions. This facilitates					
	in particular, the input of header and footer lines.					
	Program definition ( $\rightarrow$ section 5). Program position -59- is the					
	default.					
	Examples (→ section 9: Sampling operation)					

#### Checking the system using a test weight

With the built-in control of inspection, measuring and test equipment, the sequence of operations can be planned to comply with the recommendations of ISO9000, GLP (Good Laboratory Practice) and GMP (Good Manufacturing Practice). It is extremely easy to check at any time whether the LC-PVolume is operating flawlessly when connected to a balance. (First set date and time.)



Pressing the [>Conf] key normally leads to entry into the system configuration. The balance display then shows the prompt for password entry "PASS".



Skip this prompt and press the [Code] key directly. When "tare..." appears on balance display press the tare key or bar of the balance.

After taring, the following record printout starts. Load test weight and remove on completion of the printout.

	VERIFICATION
Date and time of test	13.12.94 16:25
Version of balance software	STD 10.1.19
Balance type	TYPE : AT 261
Identification number	2.00.00
Zero value	S 0.00 mg S 50000.00 mg
Value after loading test weight	5 50000:00 mg
	Ref.:
Nominal weight value, certificate	
No. or accuracy class of test weight	Act./Sign.:
Signature of test person	

# 3 General key functions/entries

Although each key of the LC-PVolume has two functions, it is not necessary to switch to a different assignment level. The system always checks the context of the key that was pressed, to decide for itself whether, e.g. a function is being called up or a number entered.



or



## Previous or next position

for selection of the program. The DeltaTrac always shows the current program position like the minute hand of a clock. In menus: Stepping up or down on the menu positions.

**Input of alphanumeric data** (for free code,  $\rightarrow$  [Code], or for character values in the configurations [like P1 or C17],  $\rightarrow$  sections 4 and 5)

[^], [v] to select A...Z/a...z, . (decimal point), - (dash) and space, [ $\downarrow$ ], or 0...9 for the next character, and [ $\downarrow$ ] to conclude the input. If the last character inputted is alphanumeric, press [ $\downarrow$ ] [ $\downarrow$ ] to conclude the input.



#### [ | key (Enter)

In (Conf) or (Prog),  $[\ \ ]$  stores the displayed value. The new entry is recorded on the printer. Otherwise the printer's paper feed advances.

The keypress is also interpreted as confirmation for various functions.



In alphanumeric entries: **Deletion of last inputted character/number,** if [] has not yet been pressed. When there is no longer a digit to the left of the cursor, the old value is again shown. Pressing the key again provides an empty input. Pressing again terminates the entry function.

The current parameter can be cleared by pressing [Clear] immediately after the parameter value is displayed. The balance display shows  $\_$ ". Confirm by pressing [ $\downarrow$ ].

Program is active:

Pressing and holding down [Clear] for a few seconds will abandon the pipette test that is currently running. Confirm 'abort' by pressing [] or refuse by pressing [Clear] again.

No program is active (standby mode):

Pressing and holding down [Clear] for a few seconds will start the procedure to delete the current program ( $\rightarrow$  section 5).



Displays the current program and its position on the DeltaTrac.

Pressing and holding down this key for a few seconds will switch from print mode to program mode.



Pressing and holding down this key for a few seconds will switch from program mode to print mode.



Starts the volumetric test using the parameters defined for the current program position.



#### [Code] key

Pressing and holding down [Code] for a few seconds while switching on LC-PVolume: The **character set** is printed out.

Entry and printout of a free **code**: The code entered must always be concluded with  $[\bot]$ .

No program is active (standby mode):

- If [Code] is pressed directly following [>Conf], a **verification** record is printed out  $(\rightarrow$  section 2).
- Pressing and holding down [Code] for a few seconds:
   A program catalog is printed out as an overview of LC-PVolume (→ section 5).

Printer mode (→section 4):
Printout of the next stable value.

# 4 Configuring the system



Pressing this key the first time enters the system configuration menu. Subsequent presses on this key after entering the correct password, switches to the next configuration parameter.

PASS

Prompt for password entry. Press  $[\ \ ]$  if no password has been defined.



Moves forward to the next configuration parameter.



Moves back to the previous configuration parameter.

Configuration parameters (characters printed in **boldface** are the current default settings):

C1	PASS	Password up to 6 digits (09, . and -). (first four characters are required for program	09
		definition and deletion; all digits for system configuration only)	

C2 La	ng Language:	1=German, 2=French,	3=English	13
-------	--------------	---------------------	-----------	----

- C4 HH.MM Time in hours and minutes, entry in 24-hour format.
- C5 nrPg Number of possible programs (test parameter 1...**60** definitions).

  (E.g. if there are only 5 programs used, set C5 to "5";

  Default position -59- is accessible only when C5 is set to "60")
- C6 ABab **O=Upper case**, 1=Mixed case letters O/1 (If set to "1", the scrolling alphanumeric character loop will also include lower case letters)
- C7 PrtM Printer mode: **0=Inactive**1=no 'TA' and no leading 'S' (stable values)
  will be printed out, 2=Print all

C8 7/8	Printer: <b>7=7 bit, even parity, no handshake,</b> 7/8 <b>baud rate=2400</b> , 8=8 bit, no parity, Xon/Xoff, baud rate=2400	
	Note: Restart LC-PVolume to activate new settings of C7/C8.	
C9 LCo	Number of line feeds after code input (3). 0255	
C10 LOt	Number of line feeds after miscellaneous 0255 printouts (3).	
C11 bUn	Barometer unit: <b>1=mmHg</b> , 2=mbar, 3=hPa 1/2/3	
C12 reJ	Printout of rejected values: 0=no printout, 1=printout of total number of rejected values, 2=printout of total number and of all (max. 30) individual rejected values.	
C13 doo	Automatic door mode (AT balance): 0=none, 0/1/2/3 1=open, 2=open/close, 3=close	
C14 eva	Working with evaporation (blank) test: 0=no, 1=yes 0/1	

To print out configuration press [Code] when the balance display shows the configuration parameter "C1 Pass".



Quits the system configuration when the title of a configuration parameter ("C..") is on the display. Otherwise, press  $[\tilde{} \tilde{} \tilde{$ 

# 5 Defining the program (test parameter)



or



Selection of the desired storage position of a new program (Display: -n- for empty positions, otherwise program name). The positions are shown by the DeltaTrac.



0...59



Goes directly to the position specified.



Displays the current program and its position on the DeltaTrac. Pressing and holding down this key for a few seconds will switch from printer mode back to standby mode.



Pressing this key the first time enters the test parameter menu. Subsequent presses on this key, after entering the correct password, switches to the next test parameter.

PASS

Prompt for password entry. Press  $[\ \ \ ]$  if no password has been defined.



or



[\lambda] Moves forward to the next test parameter (same as >Prog)

[v] Moves back to the previous test parameter

After choosing one of the test parameters (e.g. 'P1' for instrument name, [see following test parameters]), press [ $\downarrow$ ] to check the actual value. Press [ $\downarrow$ ] or [Clear] to leave it unchanged, or input your new value and press [ $\downarrow$ ] to confirm.

Test parame	ters (characters printed in <b>boldface</b> are the default settings):	Special header and footer lines (for P8P13):
P1 Inst	Name of the volumetric equipment, up to 20 characters $[0][9]$ $[\times]$ ; $[\times]$ for A to Z and space, $[\times]$ for the next character, $[\times]$ to conclude the input. Important: There has to be an input for this parameter, otherwise LC-PVolume will cease to continue.	Date/Time:  Entry: .1. Printout: 13.12.94 16:25:00 Entry: .2. Printout: 13.12.94 - 16:25 Entry: .3. Printout: 13.12.94 Entry: .4. Printout: 16:25 Entry: .5. Printout: 16:25:00
P2 Opr	Name of the operator, up to 15 characters. Valid characters are the same as in P1.	Entry: .6. Printout: Entry: .7. Printout: ====================================
P3 no.	Identification number (additional info),up to 20 characters. Valid characters are the same as in P1.	.1. Date and time (long format) .2. Date and time (short format) .3. Date
P4 noM	Nominal volume to test [ $\mu$ l]: <b>0.0</b> 0.01 10'000'000	.4. Time (short format) .5. Time (long format)
P5 -n-	Sample size: 4=Short, <b>10=Normal</b> , 4/10/30 30=Long	.6. Single line .7. Double line
P6 Cyt	Cycle time in seconds. <b>15 s</b> 199	Date/Time/Position No.:  Entry: .11. Printout: -13.12.9416:25:00- 7-
P7 Eco	EconoMode: <b>0=with instructions</b> , 1=Normal, 0/1/2 2=Compressed (see sample printouts → section 9: Sampling operation)	Entry: .12. Printout: - 13.12.94 - 16:25 7- Entry: .13. Printout: 13.12.94 7- Entry: .14. Printout: 16:25 7- Entry: .15. Printout: 16:25:00 7-
P8 †1 P9 †2 P10 †3	Header, line 1 date/time (see opposite), or 24 char. Header, line 2 date/time (see opposite), or 24 char. Header, line 3 date/time (see opposite), or 24 char.	Entry: .16. Printout: 7- Entry: .17. Printout: ======== 7-
P11 E1 P12 E2	Footer, line 1 date/time (see opposite), or 24 char. Footer, line 2 date/time (see opposite), or 24 char.	Entries .11.,, .17.: Same as entries .1.,, .7., but in addition, a position number is indicated on the right hand side of each line.

date/time (see opposite), or 24 char.

P13 E3

Footer, line 3



Quits and saves test parameter definitions.

#### Printing out program catalog



Pressing and holding down [Code]: A program catalog will be printed out as an overview of all positions.

Only the following parameters are printed out for every occupied storage position: Instrument, operator, number, nominal volume and sample size.

#### **Printing out test parameters**

For a printout of all test parameters of an instrument, select an instrument using the  $[\land]$  and  $[\lor]$  keys.



Entry into the database of the selected test. Enter password, then press  $[\ \ ]$ .



If the [Code] key is pressed while the balance display shows "P1 Inst", all **test parameters** of the selected test are printed out.

#### **Deleting test parameters**



The current instrument can be deleted from the memory by pressing and holding [Clear]. The balance display then shows "Clr.Prg". Confirmation with  $[\ \ \ \ ]$ . If a password has been defined, the balance will then ask for the password. Enter password, then press  $[\ \ \ \ ]$ .

#### Clearing individual string parameters



The current parameter can be cleared by pressing [Clear] immediately after the parameter value is displayed. The balance display shows  $\tilde{}$ \_'. Confirmation with [ $\downarrow$ ].

#### Default position: -59-

The parameter values of storage position -59-, if defined, will serve as default for all other test positions that are yet to be defined. In this case, the previous default values, as marked in this chapter are overwritten by the values of position -59-.

In the definition of test parameters for the default position, It is not necessary to enter the instrument name before moving on to the other parameters.

soon as the balance gets a stable value.

#### 6 How does LC-PVolume work?

For the blank tests, remove the beaker b1 15... Following is a step-by-step procedure on how to properly run LC-PVolume. from the balance, wait for a few seconds, ...b1 0 and return the beaker again on the Step No. Kev Display Procedure balance. Select an instrument using the  $\lceil \land \rceil$  and  $\lceil \lor \rceil$ Note: For the blank tests, no water is keys. Prepare the test instrument. added into the beaker. 2 RUN LC-PVolume is setting up for the 8 -0.31 The weighing result (mg) is displayed on Start volumetric evaluation using the test the balance. It shows the amount of parameters defined for the current storage evaporated water during the countdown position. time. 3 **tFMP** Enter the temperature reading (15... 9 **TIMER** If the current individual test is acceptable. Accept 30°C), then press [⊿]. press [Accept]. The weight result will be If a test has been conducted prior to this printed out on the test record. If an error one, the temperature entered in that test occurs during the test cycle, press was kept in memory. To view this value, [Reject]. The weight result will be Reject press [4]. Key in a new value if the discarded and the sample counter will be 6 previous temperature is not the one reset. As soon as either [Accept] or desired then press [4] again. [Reject] is pressed, LC-PVolume will again prompt the user to press [Timer]. barP Enter the barometer reading, then press Timer [괴]. Do the same procedure as in 10 Wait for the countdown display to appear. 5 temperature in case a barometer reading has been entered for a prior test. 11 -1-15... For the sample tests, remove the beaker Place a sample beaker containing water from the balance, dispense water from the on the balance. pipette into the beaker, and put the beaker back on the balance. **Timer TIMER** LC-PVolume is ready for the user to press [Timer]. WAIt The countdown display will appear as 6

Step No.	Key	Display	Procedure
12		78.13	The weighing result (mg) is displayed on the balance.
13	Accept 4 or Reject 6		Like the blank tests, each sample test could be accepted or rejected.
14			Repeat Steps 6 to 13, with a blank test being done every 10th sample test, until the last of the sample/blank tests are done.
			Note: For the short test, evaluation ends after the 4th sample test; for the normal test, after the 2nd blank test; and for the long test, after the 4th blank test.
15		tEMP	The end of the evaluation is signaled by the prompt for the second temperature reading ( $\rightarrow$ step 3).

After the second temperature is entered, LC-PVolume prints out the calculated mean volume and the results of the statistical computations (imprecision and inaccuracy).

# The Gravimetric Analysis

Before performing the test procedure, make sure that the temperatures of the equipment to be evaluated, the room, and the test liquid are identical, and as stable as possible  $(+/-0.5^{\circ}C)$  for at least 2 hours before and throughout the evaluation period.

#### Measurement of evaporation (C14=1)

This test method makes use of the weights of a series of weighing samples of a test liquid (distilled water), corrected for evaporation. Before the actual test is done, a weighing vessel (e.g. beaker) containing a small amount of test liquid is placed on the balance. Following this, the temperature and barometer pressure readings should be recorded.

In order to obtain complete and correct test sampling results, perform the following procedures: Start the timer, remove the weighing vessel, wait for a few seconds, and place the vessel back on the balance. The weighing result is the weight of the evaporated test liquid. Repeat the procedure again, but this time, add test liquid to the vessel before putting the vessel back on the balance. The weight result is the weight of the current sample.

Repeat these procedures, doing the blank test every 10 sample tests, until the last of the test has been done. Record the temperature again.

#### Preventing evaporation (C14=0)

If the test is performed using a humidity trap or is within a saturated environment, then the blank tests are no longer necessary. Setting C14 of the system configuration to 0 will omit the blank tests.

# 7 Error messages during operation of LC-PVolume

Below are error messages that may appear as a result of an incorrect operation of the program:

noProg The balance will display this error message if the program for

the current storage position is not defined.

Wrong Unit This message will be printed out if the balance is configured to a

weighing unit other than mg or g.

-RANGE- The balance will display this error message if input for

temperature or barometer readings are beyond the range for the

gravimetric analysis.

For temperature, range is from 15°C to 30°C.

For barometer pressure, range is from 600 mmHg to 800

mmHg; 800 mbar to 1067 mbar; 800 hPa to 1067 hPa.

noNom The balance will display this error message if no nominal

volume has been defined for the current program.

#### 8 What if...?

..the printer reports Entry outside the designated definition range, Input not valid? e.g. blank instrument, wrong date format, etc.

...the printer reports Wrong password or non-modifiable

Access denied? parameter.

...the system automatically No key has been pressed for at least 10

quits a function? minutes.

...the printer reports
Transmission error?

Transmission rate, number of data bits or parity is incorrectly set on the balance.

In program mode: 2400 baud, 7 bits, even parity, no handshake

2400 baud, 7 bits, even parity, no

In printer mode: handshake

(selectable in C8) 2400 baud, 8 bits, no parity, handshake

Xon/Xoff

...the printer reports
Battery discharged?

This message appears every 2 hours when the capacity of the batteries is no longer sufficient to store the data if a power outage occurs.

**Warning**: Do not switch off LC-PVolume Controller and do not disconnect from the power supply. Otherwise, the data could be lost. Replace batteries as described in Section 10: Installing the batteries.

...the printer reports
Memory reset?

The memory has been cleared due to low battery voltage. All data has been lost.

...the printer reports
Printer mode?

The LC-PVolume is in printer mode at start-up (configured in C7) or [1] was pressed and held for a few seconds. Only data from a master weighing system will be printed out by the printer. Press and hold down [Actual] for a few seconds to switch to LC-PVolume mode.

# 9 Sampling operation (examples)

System configuration: Default settings, date and time must be defined; Printout of rejected values = 0

Program definition: Nominal volume = 80  $\mu$ l; Sample size = 4; EconoMode = 0

→ section 4: Configuring the system

→ section 5: Defining the program

Keys pressed/weighings		Balance display	Printout
Program selection with keys [ $\land$ ] and [ $\lor$ ].	P 80	Position indicator with the DeltaTrac.	- 29.03.95 - 14:28 0- - THE QUALITY COMPANY Operator Dennis Lauzon
Start the volumetric evaluation with [Start].	RUN		
Enter the temperature reading. [2] [1] [.] [5] []	tEMP ⇒	21.5	Inst P 80 No. 569
Enter the barometer reading.  [7] [5] [4]	barP ⇒	755	Nominal 80.00 $\mu$ l Cycle time 15 s
Start countdown with [Timer].	tIMER	DeltaTrac indicates which sample is about to be processed.	T1 = 21.5 °C Air press. = 755.0 mmHg
Counting down. Make blank test: Remove beaker from the balance and return it back on the balance before timer	WAIt	about to be proceeda.	<pre>* Place beaker with * water on the balance. * Press <timer> when</timer></pre>
reaches 0.	b1 15 ↓		* ready for blank test.
Do not dispense water into the beaker.	b1 1		Sample # Weight
The blank test result is displayed (should be a negative	-0.31		e 1   -0.31 mg
value). Print it out on the record with [Accept].  As soon as [Accept] is pressed, the "tIMER" prompt appears again on the display.	tIMER		

# Continued...

Keys pressed/weighings		Balance display	Р	rintout
Repeat the preceding steps but instead of just returning the beaker back on the balance, dispense the contents of one pipette into the beaker for every test cycle.	78.13 77.94 77.95 77.25	Test indicator with the DeltaTrac.	1 2 3 4	78.13 mg 77.94 mg 77.95 mg 77.25 mg
After the 4 test cycles have been done, enter the temperature reading again. [2] [1] [.] [7] [.]	tEMP ⇔	21.7	T2 Mean T Mean vol	•
Automatic printout of results.			Inaccuracy E E% Imprecision	(mean error): -1.623 μl -2.03 %
			s (est.) CV(est.) Range	0.221 μl 0.28 %

System configuration: Default settings, date and time must be defined; Printout of rejected values = 1

Program definition: Nominal volume =  $50 \mu$ l; Sample size = 10; EconoMode = 1

→ section 4: Configuring the system

→ section 5: Defining the program

Keys pressed/weighings	Balance display		Prir	Printout	
Program selection with keys [\times] and [\times].  Start the volumetric evaluation with [Start].	P 50 RUN	Position indicator with the DeltaTrac.	- THE QUALIT	14:26 0- Y COMPANY Dennis Lauzon	
Enter the temperature reading.	tEMP ⇒	21.0	Inst	P 50	
[2] [1] [.] [0] [↓] Enter the barometer reading. [7] [5] [5] [↓]	barP ⇒	755	Nominal Cycle time	50.00 μl	
Start countdown with [Timer].	tIMER	DeltaTrac indicates which sample is about to be processed.	T1 = Air press. =	21.0 °C 755.0 mmHg	
Counting down. Make blank test: Remove beaker from the balance and return it back on the balance before timer reaches 0. Do not dispense water into the beaker.	WAIt b1 15 J	3.300. 10 20 p. 3000.	Sample # e 1	Weight -0.22 mg	
reaches of Do not dispense water into the beaker.	b1 1		1 2	53.07 mg 52.61 mg	
The blank test result is displayed. Print it out on the record with [Accept].	-0.22		3 4 5	50.98 mg 51.88 mg 51.16 mg	
	шмер		6	51.16 mg 51.97 mg 51.61 mg	
As soon as [Accept] is pressed, the "tIMER" prompt appears again on the display.	tIMER		8	51.95 mg	
			9 10	50.58 mg 52.06 mg	

# Continued...

Keys pressed/weighings		Balance display		Printout
Repeat the preceding steps but instead of just returning the beaker back on the balance, dispense the contents of one	53.07 52.61	Test indicator with the DeltaTrac.	e 2	-0.51 mg
pipette into the beaker for every test cycle.	50.98 51.88		T2 Mean T	= 20.8 °C 20.9 °C
To redo test cycles with undesirable results, press [Reject] rather than [Accept] when the weight result is displayed.	51.16 51.97 51.61		Mean vol	
[Rejected value]	0.00 51.95		E	y (mean error): $2.311 \mu l$
[Rejected value]	100.23 50.58		E% Imprecision	4.62 %
[Rejected value]	75.23 52.06		s CV	0.748 $\mu$ l 1.43 %
After the 10 test cycles have been done, perform another blank test.	-0.51		Range Rejected	2.50 μl values: 3
After the second blank test, enter the temperature reading again. [2] [0] [.] [8]	tEMP ⇒	20.8		14:28
Automatic printout of results, including total number of rejected values.				

System configuration: Default settings, date and time must be defined; Printout of rejected values =2

Program definition: Nominal volume =  $100 \mu l$ ; Sample size = 30; EconoMode = 2

→ section 4: Configuring the system

→ section 5: Defining the program

Keys pressed/weighings		Balance display	Pr	intout
Program selection with keys [^] and [v].  Start the volumetric evaluation with [Start].	P 100 RUN	Position indicator with the DeltaTrac.	- THE QUALI	- 14:20 0- TY COMPANY Dennis Lauzon
Enter the temperature reading.  [2] [0] [.] [7] [  Enter the barometer reading.	tEMP ⇒ barP ⇒	20.7 758	Inst No. Nominal	•
Enter the barometer reading. [7] [5] [8] [↓]	buir 🛶	700	Cycle time T1 Air press.	= 20.7 °C
Start countdown with [Timer].  Counting down. Make blank test: Remove beaker from the balance and return it back on the balance before timer reaches 0.	tIMER  WAIT b1 15  U b1 1	DeltaTrac indicates which sample is about to be processed.	e 1  1 2 3 4 5 6 7	-0.28 mg 98.18 mg 98.25 mg 99.70 mg 95.88 mg 99.98 mg 96.18 mg 98.22 mg
Do not dispense water into the beaker.  The blank test result is displayed. Print it out on the record with [Accept].	-0.28		8 9 10 e 2 11 12 13	99.29 mg 99.35 mg 100.07 mg -0.53 mg 99.17 mg 99.24 mg 99.69 mg


# Continued...


Keys pressed/weighings	Balance display		Printout
As soon as [Accept] is pressed, the "tIMER" prompt appears again on the display.	tIMER	14 15 16	99.65 mg 99.88 mg 102.00 mg
Repeat the preceding steps but instead of just returning the beaker back on the balance, dispense the contents of one pipette into the beaker for every test cycle.	98.18 Test indicator with the 98.25	17	99.95 mg 99.63 mg 101.58 mg 99.21 mg
After the 10 test cycles have been done, perform another blank test.	-0.53	e 3 21 22 23	-0.49 mg 100.19 mg 98.35 mg 100.42 mg
Perform the cycle tests for the next 10 test cycles. Press [Reject] to redo the tests with undesirable results.	99.17 Test indicator with th 99.24 99.69	26 27	99.39 mg 100.11 mg 100.53 mg 100.83 mg
[Rejected value]	97.46 99.65 99.88	28 29 30 e 4	100.36 mg 100.27 mg 99.93 mg -0.40 mg
[Rejected value]	89.75 102.00 99.95	T2 Mean T	= 20.9 °C 20.8 °C
	99.63 101.58 99.21	Mean vol Inaccuracy	100.243 μl y (mean error):
Perform another blank test.	-0.49	E E%	0.243 μl 0.24 %


# Continued...

		Balance display	FII	ntout
Perform the last 10 test cycles. Press [Reject] to redo the tests with undesirable results.	100.19 98.35 100.42 99.39	Test indicator with the DeltaTrac.	Imprecision: s CV Range	1.295 μl 1.29 % 6.14 μl
[Rejected value]	-0.22 100.11 100.53 100.83 100.36 100.27 99.93		Rejected val	97.46 mg 89.75 mg -0.22 mg 99.64 mg
After the last 10 test cycles, the last blank test is done.	99.64			
An error was made during the last blank test. Press [Reject] and redo the test.	-0.40			
Enter temperature reading again. [2] [0] [.] [9] [↓]  Automatic printout of results, including rejected values.	tEMP ⇔	20.9		

## 10 System startup and maintenance

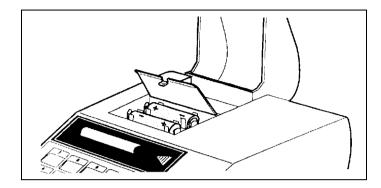

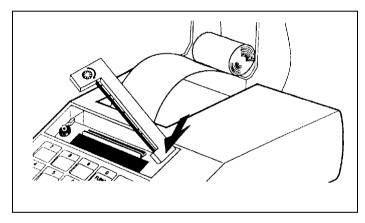

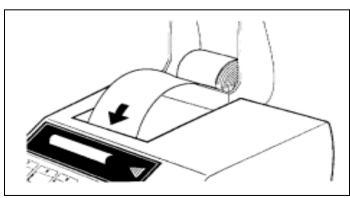

#### Installing the batteries

The two enclosed batteries power the clock and the data memories for the settings, register data and statistics when the LC-PVolume Controller is not connected to the power supply.

**Please note:** If work has already been performed with the system and the existing data have to be retained, the LC-PVolume Controller must be connected to the power supply and switched on when the batteries are changed. Otherwise, the data will be lost within a few minutes.

- Raise paper cover and pull paper backward out of printer
- Open battery compartment and insert batteries

#### Replacing the ribbon


- Remove paper from printer
- Remove ribbon cover
- Remove ribbon cartridge
- Insert new ribbon cartridge (tighten with the tensioning wheel if necessary) and reattach ribbon cover

#### **Inserting paper**

- Insert paper in the slot of the printer in the direction of the arrow
- Switch on LC-PVolume Controller (toggle switch at rear right)
- Press and hold the 「↑] key until sufficient paper has been drawn in

#### **Connecting LC-PVolume to balance**


Connect LC-PVolume to the balance using the appropriate cable (→ Appendix B). Use one of the two small, round and green CAN plugs to connect a CAN balance, or the 9-pin RS232 plug for all other balances.

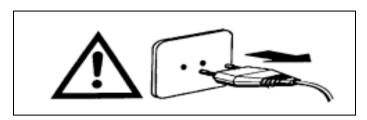

#### **Cautionary note:**


If the printer or the power fuse has to be replaced, the housing must be opened.

Before opening the instrument, make sure that the power cable is unplugged!

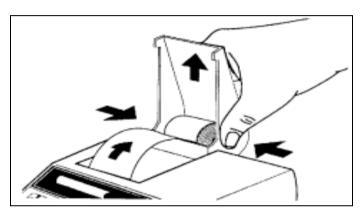

#### Replacing the printing unit

The printing unit can be replaced. Order number in Appendix C: Standard and optional equipment. The instructions for replacing the printing unit are enclosed with the printer.

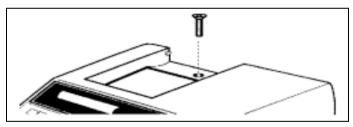


#### Replacing the power fuse

- Pull paper out of printer and remove roll
- Gently press paper cover together until it disengages from the controller housing
- Remove paper cover
- Unscrew housing screw
- Carefully open housing (pay attention to ribbon cable connection to top housing; the screening plate has been omitted from the illustration)

Replace fuse
 230V: slow-blow 100mA, 250V
 115V: slow-blow 200mA, 250V

• Close housing and tighten screw

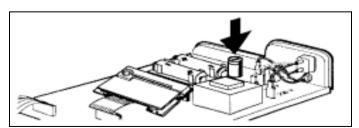

### Appendix A: Technical data of LC-PVolume Controller

#### **Printing unit**

Type of printing unit Dot-matrix printer 5x9 dots, 24-character/line

Epson/IBM character table No. 4

Printing speed Faster than 1 line per second

Ribbon cartridge Exchangeable, black

Paper roll Normal paper 58 x  $\varnothing$  51 mm, integrated in

housing, commercial size

**Clock** Time and date, EURO or US format representation,

leap year taken into account

Interface RS232C bi-directional, 2400 baud fixed

LC-PVolume mode: printer mode selectable:

Number of data bits77or8ParityevenevennoneHandshakenonenoneXon/Xoff

Printer buffer 128 characters

Mechanical port

Temp. range/humidity 0...40°C/15...85% r.h.

Approvals EN60950 (equivalent to UL1950), CSA

**Power supply** 

Line voltage, 100...120V, 160mA or 220...240V,

Consumption, frequency 80mA, 50...60Hz

**Dimensions,** W x D x H = 157x210x85mm, weight net 1.5kg (incl. paper roll)

# Appendix B: Standard and optional equipment

# Included in the standard equipment of the METTLER TOLEDO LC-PVolume:

[One of them is included]

Order No.

1 Paper roll	→ below
1 Ribbon cartridge, black	→ below
2 Batteries UM3 (1.5V) accord. to IEC LR 6	commercial
Fuse 230V (Slow-blow 100mA, 250V) or	_
Fuse 115V (Slow-blow 200mA, 250V)	
according to IEC 127 / III	commercial
Operating Instructions German	21900418
Operating Instructions French	21900419
Operating Instructions English	21900302

#### **Optional accessories:**

Paper rolls, set of 5 Ribbon cartridges, black, set of 2	72456 65975			
Connection cable for AE (Option 012)	59759			
Connection cable for LC-PVolume - Bo (AT, MT, UMT, AM, AJ Option 018)	229029			
Connection cable SQC14 – balance RS232, (PG-S, PB, Spider1)		11101051		
For CAN/SICS balances: Connection/extension (max 10 m) cable				
for LocalCAN, 0.3 m for LocalCAN, 1.0 m	LC-LC03 LC-LC1	239270 229161		

# Index

A	G	Power fuse21 Previous position5, 6
Aborting5	Gravimetric Analysis3, 11	Printer mode
Accept key10, 14		Prog. key9
Actual key5, 7		Program catalog
Air pressure10, 13	Humidity trap11	R
Alphanumeric keys6	I	-
В	Identification number8	Record footer8 Record header8
Barometer unit7	Imprecision11	Reject key10
Batteries20	Inaccuracy11	Rejected values
Blank test10, 13	Input of alphanumeric code5	Ribbon20
C	L	S
Character set6	Language6	Sample size8
Clear5, 9	LC-PVolume Controller22	Saturated environment11
Code key4, 6, 9	Line feeds7	Special header and footer lines8
Configuration4	M	Start key10, 13
Configuration parameters6	Mean volume11	System configuration6, 7
Cycle time8	would volume	T
D	N	Temp. reading10, 11, 13, 14, 19
Date and time4, 6, 8	Name of operator8	Test parameters
Door (mode for AT balance)7	Name of volumetric equipment8	Timer key
F	Next position	V
	Nominal volume8	V
Econo Mode8	P	Verification4, 6
Enter key5 Error messages12	Paper insertion on printer20	
Evaporation test (active/inactive)	Password6	

To protect your METTLER TOLEDO product's future:
METTLER TOLEDO service assures the quality, measuring accuracy and preservation of value of all METTLER TOLEDO products for years to come.
Please send for details about our attractive terms of service.
Thank you.



© Mettler-Toledo GmbH 1995

ME-21900302A Printed in Switzerland 9911/2.12

**Mettler-Toledo GmbH,** CH-8606 Greifensee, Switzerland, Phone +41-1- 944 22 11, Fax +41-1-944 34 30, Internet: http://www.mt.com Subject to technical changes and to the availability of the accessories supplied with the instruments.