
METTLER TOLEDO

Emulation B → M for B series balances **Emulation B → M für Waagen der B-Linie** **Emulation B → M pour les balances des ligne B**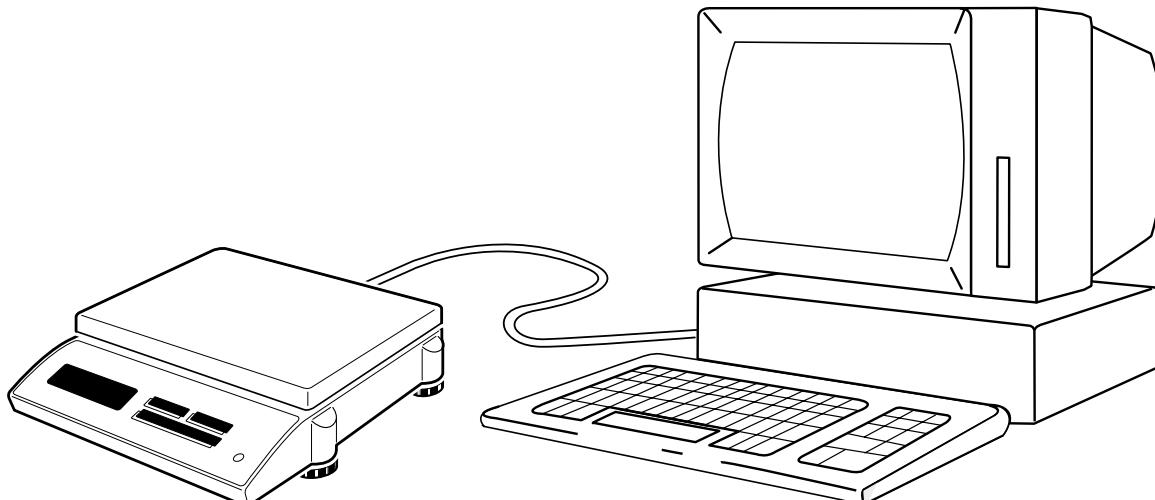

English

Introduction

The „Emulation B → M“ makes the B balance line (PB, SB) compatible with the M and J series. „B → M“ emulates the menu, functions and the interface of the M and J balances on the B balances. This emulation facilitates integration of the B balances in existing systems which interchange data with a host computer or which are controlled by a computer as the existing programs do not have to be modified. The „B → M“ software is accommodated on an EPROM.

Standard equipment

- EPROM „Emulation B → M“ (with RS232C interface: EPROM for balance board; with LocalCAN interface: OTPROM for LC-B option board).
- Operating instructions “DataPac-M” (in English only)
- Installation information “Emulation R/G → M”
- Installation “Emulation B → M” (this document)

Installation and settings

- For balances with a LocalCAN interface, you need the set 21301729. Replace the OTPROM on the option board by the OTPROM supplied.
- For balances with an RS232C interface, you need the set 21301730. Replace the EPROM on the balance board by the EPROM supplied.

For the setting of the interface parameters, please consult the installation information „Emulation R/G → M“. For the menu settings, see page 18: The „Emulation R/G → M“ on G-S balances.

Important: With the B balance, the CAL function appears before entry into the menu.

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