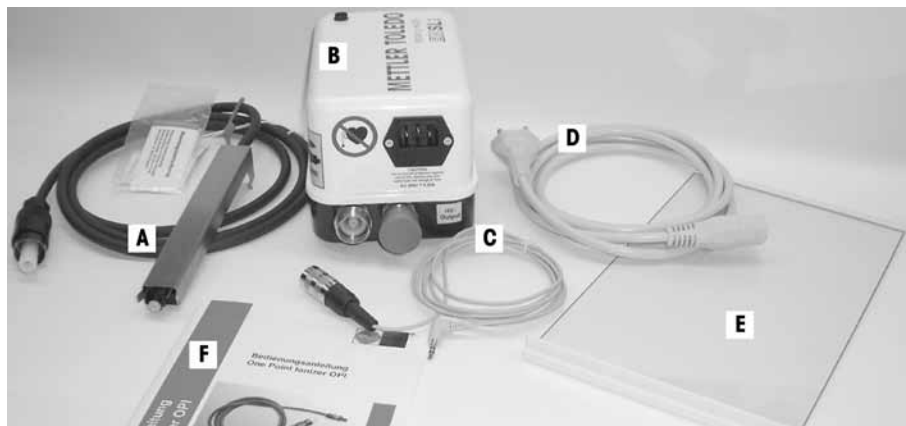
Installation Instructions **Antistatic Kit for XP Analytical Balances**
 Montageanleitung **Antistatik Kit für XP Analysenwaagen**
 Notice de montage **Kit antistatique pour les balances d'analyse XP**
 Instrucciones de montaje **Kit antiestático para balanzas analíticas XP**
 Istruzioni di montaggio **Kit antistatico per bilance analitiche XP**

Components

- | | |
|-------------------------------|---------------------------------|
| A Single-point ionizer | B AC adapter |
| C Control cable | D Power supply cable |
| E Side door | F Operating Instructions |

Note

Position the single-point ionizer on the left or right side of the balance.
The AC adapter can have two single-point ionizers connected to it.

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Italiano

**CAUTION****Damage of Device**

The power cable used to connect the ionizer must have a ground wire and be plugged into an outlet with a functioning ground connection.

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**DANGER****Persons with cardiac pacemakers**

Persons with cardiac pacemakers must always maintain a safety distance **of more than 50 cm** from the ionizer.

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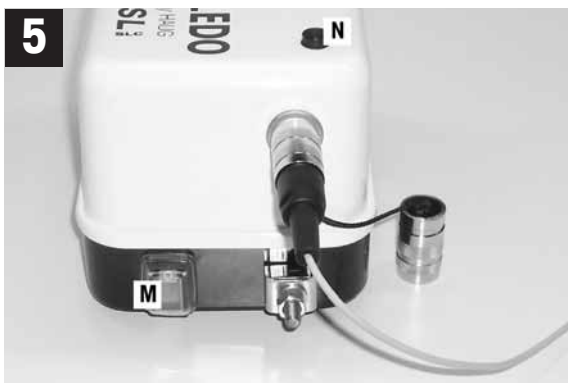
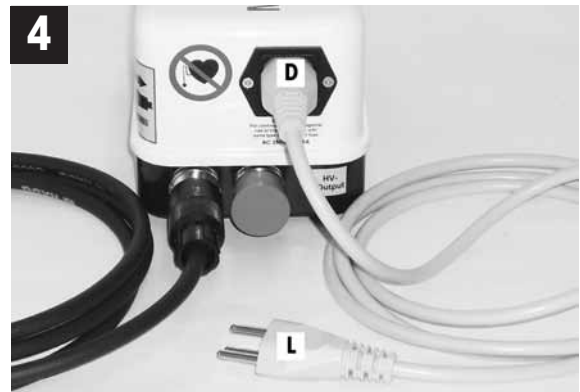
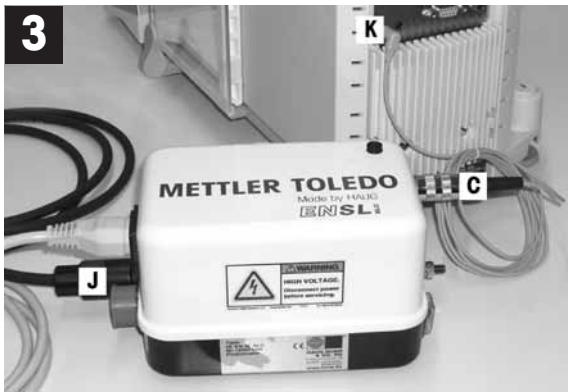
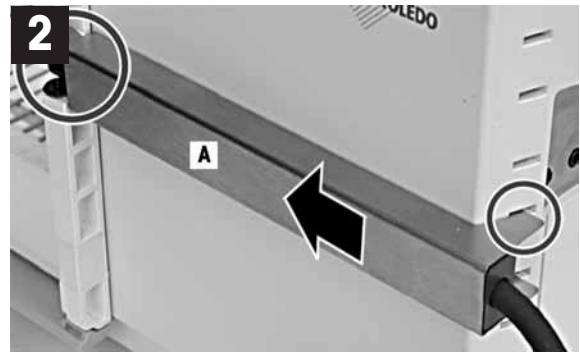
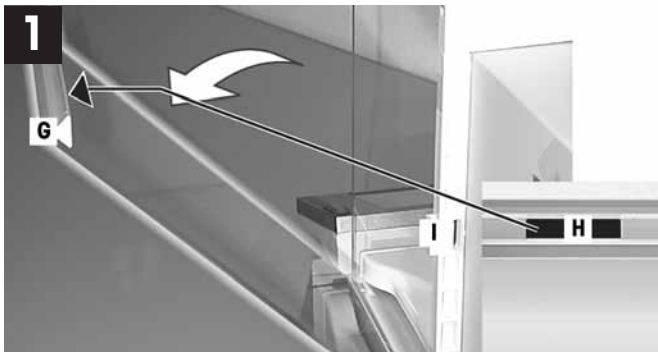
English

- 1 • Swap the side door (G) for the door (E) marked with "Conducting Glass" (H) supplied with the antistatic kit.
 - Pull the middle clip (I) (recommended position) off.
- 2 • Push the point ionizer (A) in from the back. If necessary, slightly bend the strips so that the ionizer is held firmly.
- 3 • Screw the ionizer cable (J) tightly onto the AC adapter.
 - Screw in the control cable (C).
 - Plug the control cable into Aux 1 or Aux 2 (K).
- 4 • Plug in power cable (D).
 - Connect power cable (L) to power supply.
- 5 • Switch On/Off (M).
 - ⇒ Control lamp (N) lit when the ionizer is active.

Information about settings for the antistatic kit is contained in the operating instructions for XP analytical balances.

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© Mettler-Toledo AG 02/2012
Mettler-Toledo AG