English

## Installation instructions

BT-BLD Auxiliary display, 11132555

Deutsch

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Français

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Afficheur auxiliaire BT-BLD, 11132555

Español

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Display auxiliar BT-BLD, 11132555

Italiano

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Indicatore ausiliario BT-BLD, 11132555

METTLER TOLEDO

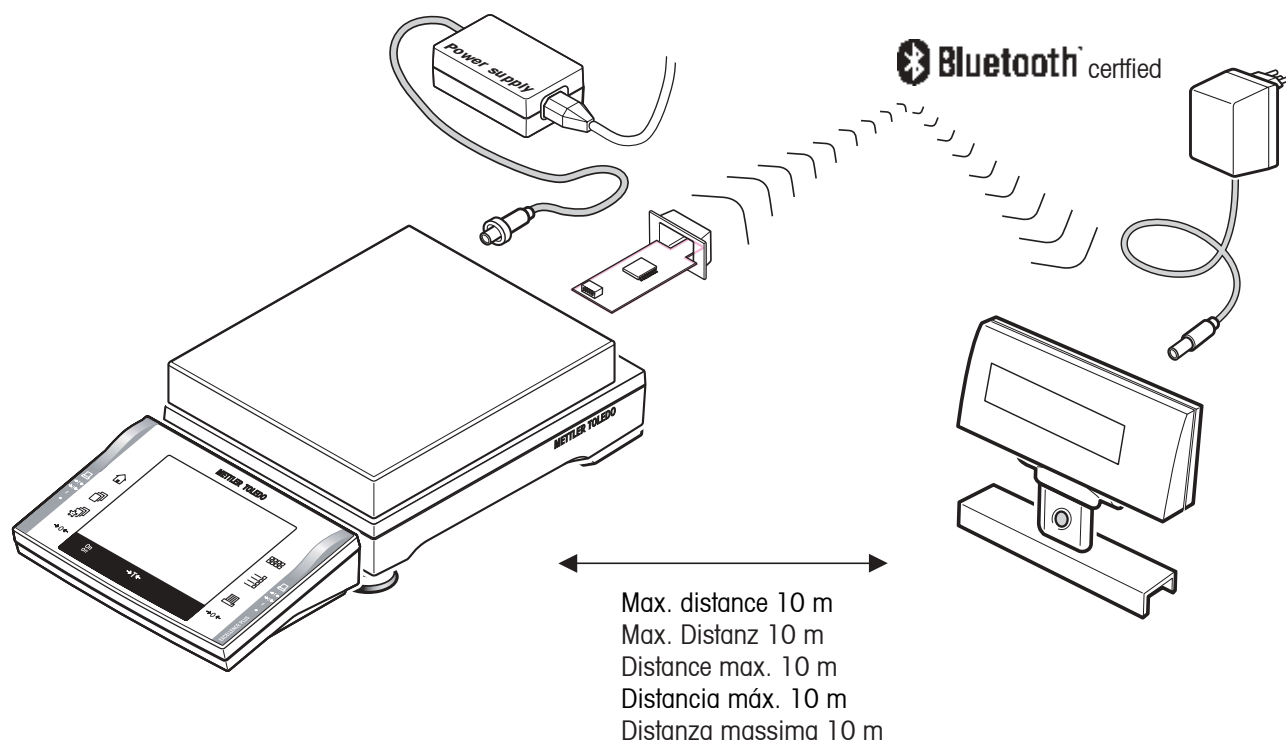
BT-BLD auxiliary display for wireless connection to an XP/XS precision balance.  
(BT or BTS interface option required)

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
Bluetooth is a registered trademark of Bluetooth SIG, Inc.

© Mettler-Toledo GmbH 2004 11780629A Printed in Switzerland 0409/33.45

Mettler-Toledo GmbH, Laboratory & Weighing Technologies, CH-8606 Greifensee, Switzerland  
Phone +41-44-944 22 11, Fax +41-44-944 30 60, Internet: <http://www.mt.com>

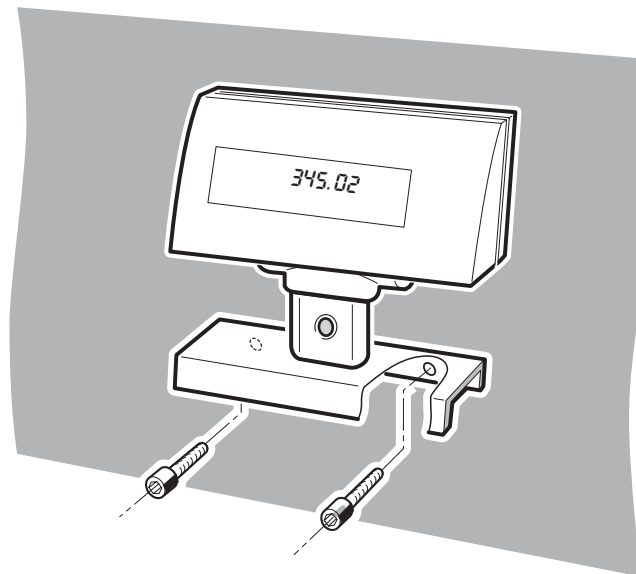Wall mounting  
Wandmontage  
Montage mural  
Montaje en pared  
Montaggio a parete

#### Safety warning

- If equipped with the power supply unit delivered ex works, balances fitted with the BT-BLD auxiliary display must **not be operated in potentially explosive environments**, and only indoors in a dry environment.
- Before connecting the power supply unit, check that the voltage marked on it is identical with the voltage of the local power supply. If this is not the case, please contact your local MT dealer.
- Use only a tested AC adapter with SELV output current.

#### Connecting the auxiliary display to the balance

- Place the auxiliary display within receiving distance to the balance. Maximum distance: 10 m
- Connect the device in accordance with the Installation and Operating instructions for the BT or BTS interface option. As soon as the device to be connected has been set up for the BT-BLD, the auxiliary display is ready to operate. The connection may take 10 to 20 seconds to establish.

The connection status is indicated by the blue Bluetooth lamp (1):

- Lamp flashing: No connection
- Lamp continuously lit up: Connection present

#### Technical data

Standard equipment: Power supply to auxiliary display, Input 12–16 VAC, 50/60 Hz, 3 VA or 9–22 VDC, 3 W

Country-specific power unit:

Output 12 VAC, 500 mA

– Euro	230V/50Hz/80mA	11103740
– Euro/Schuko	230V/50Hz/80mA	11103744
– UK	240V/50Hz/80mA	11103742
– USA	120V/60Hz/10W	11103741
– Japan	100V/50Hz/10W	11103743

Output 12 VAC, 1.0 A

– Universal	220–240V/50Hz/100mA	11103745
-------------	---------------------	----------

(Bench version, requires additional country-specific cable)

Dimensions: (W/D/H) 182/60/168 mm

Net weight: 620 g (1550 g complete with packing)

Environmental conditions: Ambient temperature 10 °C ... 30 °C  
Relative atmospheric humidity 15 % ... 80 % non-condensing  
Height above sea level ≤ 4000 m

Degree of protection: Pollution degree 2  
Overvoltage category Class II

Data transmission: Wireless Bluetooth® technology Max. range of 10 m

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