### **Operating Instructions**


# METTLER TOLEDO PL-L Line of Classic Light Balances






#### Operating instructions in a nutshell



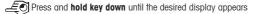




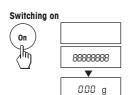


#### Switching off





#### Simple weighing

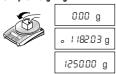


#### Adjusting (calibration) external

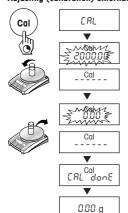

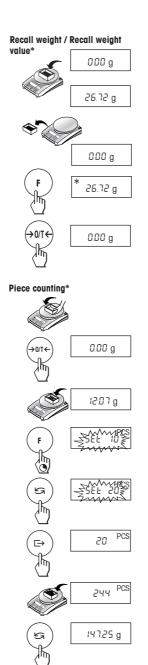

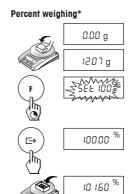

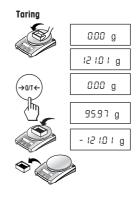





\* These functions must be activated in the menu (section 4.3.2)

#### Contents

1	Getting to know PL-L balances line	4
1.1	General	4
1.2	Layout of balances	4
1.3	Overview of key functions	5
2	Startup	6
2.1	Unpacking / standard equipment	6
2.2	Cautionary notes	6
2.3	Setting up, leveling, preparations for weighing below the balance	6
2.4	Adjusting (calibration)	8
3	Weighing	9
<b>3</b>	Weighing	
•	• •	9
3.1	On/Off switching	9
3.1 3.2	On/Off switching	9 9 9
3.1 3.2 3.3	On/Off switching	9 9 9
3.1 3.2 3.3 3.4	On/Off switching	9 9 9
3.1 3.2 3.3 3.4	On/Off switching	9 9 9 9

5	Functions	15
5.1	Recall weight / Recall weight value	15
5.2	Piece counting	16
5.3	Percent weighing	17
5.4	Switching weight units	17
6	Technical data, options,	
	optional equipment	18
6.1	Technical data	18
6.2	RS232C interface	20
6.3	MT-SICS Interface commands and functions	20
6.4	Optional equipment	22
6.5	Dimensional drawing (in mm)	23
7	Appendix	24
7.1	Typical printouts from METTLER TOLEDO	
	RS-P26 and LC-P45 printers	24
7.2	What if?	25
7.3	Maintenance and cleaning	26
7.4	Declaration of conformity	27

#### 1 Getting to know PL-L balances line


#### 1.1 General



#### **Balance features**

- The PL-L balance line comprises portable balances with a readability of 0.01 g to 1 g. The weighing ranges extend from 210 g to 6.1 kg.
- In addition to basic operations such as weighing, taring and adjusting (calibration) miscellaneous functions such as "Recall weight", "Piece counting weighing" or "Percent weighing" can be activated.

#### 1.2 Layout of balances

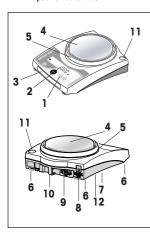

- I Keys
- 2 Display
- **3** Model plate with the following data:

"Max": maximum capacity
"d": readability

"Min": minimum capacity (recommended minimum load, only relevant for

certified balances)


"e": verification scale interval (smallest display increment tested during certification, only relevant for certified balances)

4 Weighing pan

- 5 Draft shield element (not all models)
- 6 Leveling feet (not all models)
- 7 Hanger opening for weighing below the balance (underside of balance)
- 8 AC adapter socket
- 9 RS232C interface (optional)
- 10 Lug for optional antitheft device
- 11 Leveling control (not all models)
- 12 Compartment for batteries

#### 1.3 Overview of key functions

The balances have two operator control levels: the **weighing mode** and the **menu**. The function of each individual key depends on the operator control level and how long the key is pressed.

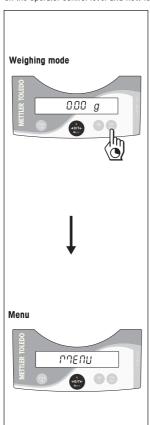

Key function	Key functions in weighing mode						
Press brie	fly 🖑	Press and hold down					
1/10d	Reduce readability	Cal • Adjust (calibrate)					
On →0/T← C	<ul><li>Switch on</li><li>Zero/tare</li><li>Cancel function</li></ul>	Off • Switch off					
S	Switch     Change settings	F • Call function; A function must be activated in the menu, otherwise "F nonE" appears in the display					
⊖	Transfer weighing data via interface with activated printer  Confirm settings	Menu • Show menu (hold key down until MENU appears)					

Key functions in menu mode					
Press briefly	, <u>(</u> m)	Press and hold down			
1/10d	Change settings Reduce value by 1 step	1/10d • Reduce value rapidly			
С	Close menu (without saving changes)	-			
S	Change settings Increase value by 1 step	• Increase value rapidly			
□ •	Select next menu item	Menu • Save changes and close menu			

#### 2 Startup

#### 2.1 Unpacking / standard equipment



The standard equipment for every balance comprises:


- · AC adapter, to national standard
- Weighing pan, Weighing pan support, draft shield element (depending on model)
- Operating Instructions
- Protective cover (placed on the balance over the weighing cell cone) with instruction sheet. This protective cover must not mislaid. It will be needed again later to protect the cone when changing batteries (underside of balance).
- In-use covers are available as optional extras (Section 6.4). In the case of models having the large weighing pan (ø 160 mm) the antistatic plate a) (secured by two screws) and the adapter ring b) must also be removed in order to fit the in-use cover.

#### 2.2 Cautionary notes



- PL-L balances must not be operated in hazardous areas with the standardsupply AC adapter.
- Before connecting the AC adapter, verify that the voltage printed on it corresponds to the local AC power supply voltage. If this is not the case, please contact your local METTLER TOLEDO dealer.
- PL-L balances may only be used indoors in a dry environment.
- For use with CSA Certified (or equivalent approved) power source, which must have a limited and SELV circuit output.

#### 2.3 Setting up, leveling, preparations for weighing below the balance

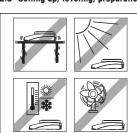


#### The optimum location

The correct location makes an important contribution to the accuracy of the weighing results of high-resolution analytical and precision balances.

- Stable, vibration-free position as horizontal as possible
- No direct sunlight
- No excessive temperature fluctuations
- No drafts

The best location is on a stable bench in a corner protected against drafts, as far away as possible from doors, windows, radiators or the louvers of air conditioners.

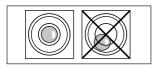

#### Levelina

Some models are equipped with a level glass and two or four leveling feet to compensate for minor irregularities in the surface on which the balance stands. The balance is exactly horizontal when the air bubble is in the middle of the level glass.

**Note:** The balance should be leveled each time it is moved to a new location.


#### Preparations for weighing below the balance

To carry out weighing operations below the balance, get rid of the special cover on the underside of the balance. (Note: never put the balance without the protective cover over its cone down on its head, only on its side!). This exposes the opening for the hanger, making weighing below the balance possible.



#### Antitheft device

All models are provided with a lug for attaching an antitheft device (see optional equipment in Section 6.4).

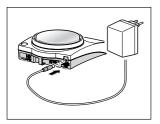

#### Power supply

- → Plug the AC adapter into the AC adapter socket on the balance, and connect to the power supply.
- → The balance performs a self-test. This test is finished when "OFF" appears.
- → Press the «On» key briefly: the balance is in operational readiness. Before any work is performed with the balance, it must be adjusted (Section 2.4).

#### **Battery operation**

Models in the PL-L line of balances can also be operated independently of the AC power supply by using their batteries. To do this, always fit the protective cover over the weighing cell cone first, then open the cover of the battery compartment on the underside of the balance and insert the batteries.

Caution: ensure correct polarity (as specified inside the battery compartment).

Close battery compartment again.

When the balance is operating on its batteries, the border around the battery symbol in the display lights up. The number of segments that are lit is an indicator of battery condition (3 = fully charged, 0 = discharged). When the batteries are almost completely discharged, the last segment flashes.

Recommended battery type: AA (LR6) 1.5 V alkali-manganese.

NiMH (nickel-metal hydride) rechargeable batteries, which are recharged in an external battery charger, can be also be used. The intervals between recharging are not as long as the service life of a nonrechargeable battery.

#### Notes

- Batteries are not included in the standard supply.
- Battery operation is automatically overridden when the AC adapter is connected to the AC power supply.
- To prolong battery (disposable or rechargeable) life, it is advisable to activate «Auto shut» in the menu (see Section 4.3.7).
- All discharged batteries must be disposed of in an environmentally responsible manner. No attempt must be made to incinerate or disassemble them.




#### 2.4 Adjusting (calibration)

To obtain accurate weighing results, the balance must be adjusted to match the gravitational acceleration at its location.

#### Adjusting is necessary

- before the balance is used for the first time
- at regular intervals during weighing service
- after a change of location

To obtain accurate results, the balance must be left switched on for 30 minutes to reach operating temperature before starting the adjustment procedure.

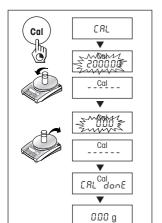

#### Adjusting with external weight

- → Have required adjusting weight ready.
- → Unload weighing pan.
- → Press and hold the «Cal» key down until "CAL" appears in the display. Release key.

The required adjustment weight value flashes in the display.

- → Place adjustment weight in centre of pan. The balance adjusts itself auto-matically.
- → When "0.00 g" flashes, remove adjustment weight.


The adjusting is finished when the message "CAL done" appears briefly in the display, followed by "0.00 g". The balance is again in weighing mode and ready for operation.

#### Notes

- Certified PL-L models cannot be adjusted by the user, because of weights and measures legislation.
- This adjustment procedure can be terminated at any time with the «C» ("Cancel") key. The balance reverts to weighing mode.

#### 3 Weighing

#### 3.1 On/Off switching

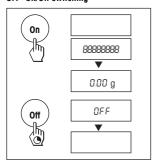

#### Switching on

→ Remove any load from weighing pan and press «On» key briefly.

The balance performs a display test (all segments in the display light up briefly).

When zero is displayed, the balance is ready for operation.

#### Switching off

→ Press and hold the «Off» key down until "OFF" appears in the display. Release the key.


#### 3.2 Simple weighing



- → Place weighing sample on the weighing pan.
- → Wait until the stability detector "o" disappears.
- → Read the result


#### 3.3 Faster weighing with reduced readability

The balance has the facility for speeding up the weighing operation by reducing its readablity (number of decimal places):

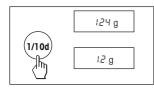

- → The balance is operating with its normal readability and speed.
- → Press the «1/10d» key and ...
- → ... the balance operates with reduced readability (one decimal place less), but displays the weighing result quicker. Pressing the «1/10d» key briefly again toggles the balance back to its full readability.

#### 3.4 Taring

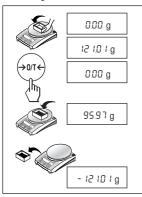

- → Place empty container on the balance.
- → The weight is displayed.
- $\rightarrow$  Press the « $\rightarrow$ 0/T $\leftarrow$ » key briefly.
- → Add weighing sample to container. The net weight is now displayed.

If the container is removed from the balance, the tare weight will be shown as a negative value.


The tare weight remains stored until the «  $\rightarrow$  0/T  $\leftarrow$  » key is pressed again or the balance is switched off.

Overview of menu

#### 4 Menu

#### 4.1 Overview

In the menu you can change the weighing unit, select additional functions and carry out various settings. A description of the individual menu options is given in Section 4.3.

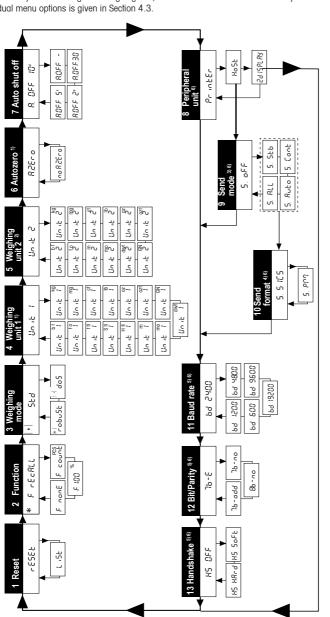

With certified balances, this menu option has a fixed setting and cannot be changed. With certified balances, only those weighing units allowed by the appropriate national

Notes

weights and measures legislation may be selected.

This menu option is only shown if "Host" has been selected in menu option 8 (Peripheral unit).

Menu option Factory setting

> These menu options are only shown if "Host" or "Printer" has been selected in menu option 8 (Peripheral unit). This menu option is only shown if "S.oFF" has not been selected in menu option 9 (Send mode). 4 2


6) Only displayed if an interface has been installed.

#### 4.2 Menu operation




#### Opening the menu

In weighing mode, press and hold down the **«Menu»** key until "MENU" appears in the display. Release the key: the 1st menu option is displayed.

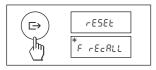

#### Select menu options

The « > » key is used to select individual menu options with their current settings one after the other.

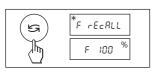

#### Change settings

Pressing the «\$\sigma\) wey displays the next setting; pressing the «1/10d» key displays the previous one. Once the desired setting appears in the display, the next menu option can be selected («\$\sigma\) or you can close the menu (see following Section).



#### Saving settings and closing the menu

Hold the **«Menu»** key down until "StorEd" appears in the display. Release the key and the balance reverts to weighing mode. All changes are saved.



#### Abort

Press the  ${}^{\rm c}$  key briefly. The balance reverts to weighing mode. Changes are  ${\it not}$  saved.

#### Note

If no entry is made within 45 seconds, the balance reverts to weighing mode. Changes are **not** saved.


#### 4.3 Description of menu options

#### 4.3.1 Reset or recording of balance settings (1st menu option "RESET")



#### Reset balance settings

→ Select "Reset", press and hold down the «Menu» key until the message "r donE" confirms that all menu settings have been reset. The balance then reverts to weighing mode and works with the factory settings (Section 4.1).

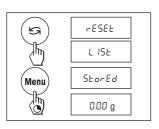


#### Recording balance settings

→ Select "List" and hold down the «Menu» key until the message "StorEd" is displayed.

The current balance settings are transmitted to the peripheral device connected to the optional RS232C interface. To do this the setting "Printer" must always be selected at the 8th menu option (Peripheral unit). The current balance settings are saved at the same time.

#### 4.3.2 Functions (2nd menu option / see Section 5 for their use)

In addition to simple weighing, the following functions can be selected with the «S» key:

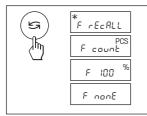

F rECALL Recall weighing
F count Piece counting
F 100 % Percent weighing
F nonE No function, simple weighing


#### 4.3.3 Weighing mode (3rd menu option)



This setting allows you to adapt the balance to the weighing mode. Select "Std" (standard) for all normal weighing processes. With "doS" (dosing) - for dispensing substances in liquid or powder form - the balance reacts very rapidly to the slightest changes of weight. With "robuSt" (absolute weighing) the balance only reacts to more significant changes in weight, so that the weighing result is very stable.

#### 4.3.4 Weighing unit 1 (4th menu option "UNIT 1")

Depending on requirements, the balance can operate with the following units (possible with certified balances only if permitted by national weights and measures legislation):

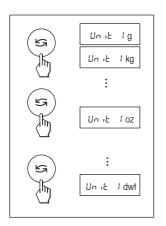

Unit	Conve	rsion factor	Comments
g gram			factory setting
kg kilogram	1 kg	= 1000 g	not with 0.1 mg and
mg milligram	1 mg	= 0,001 g	1 mg balances with 0.1 mg and 1 mg balances
ct carat	1 ct	= 0.2 g	
lb pound	1 lb	= 453,59237 g	not with 0.1 mg
			balances
oz ounce	1 oz	= 28,349523125 g	
ozt troy ounce	1 ozt	= 31,1034768 g	
GN grain	1 GN	= 0,06479891 g	not with 1 g balances
dwt pennyweight	1 dwt	= 1,55517384 g	
mo momme	1 mom	ı = 3,75 g	
m Mesghal	1 msg	≈ 4,6083 g	
H tl Hong Kong tael	1 tlh	= 37,429 g	
Stl Singapore tael	1 tls	≈ 37,7993641666667 g	The Malaysian tael has the same value
t tl Taiwan tael	1 † #	= 37.5 g	
t o tola	1 tola	= 11,6638038 g	

#### 4.3.5 Weighing unit 2 (5th menu option "UNIT 2")

If it is required to show the weighing result in weighing mode in an additional unit by pressing the « 🖘 » key, the desired second weighing unit can be selected in this menu option. The same weighing units are available as under "UNIT 1", with the exception of the toel units ("H tl", "S tl" and "t tl").

bt baht

1 baht = 15,16 g

#### 4.3.6 Autozero (6th menu option / see overview and notes in Section 4.1)

This menu option allows you to switch the automatic zero correction on or off.



#### Autozero switched on


The zero point is automatically corrected (e.g. if drift occurs or the weighing pan becomes dirty).

#### Autozero switched off

The zero point is **not** automatically corrected. This setting is advantageous for special applications (e.g. evaporation measurements).

#### 4.3.7 Auto shut off (7th menu option)

If the automatic shut off function is activated, the balance automatically switches itself off after a selected period of inactivity (i.e. with no key being pressed or changes of weight occurring):

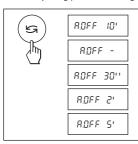

A.OFF 10' Automatic shutoff after 10 minutes inactivity
A.OFF - Automatic shutoff **not** activated
A.OFF 30' Automatic shutoff after 30 seconds inactivity
A.OFF 2' Automatic shutoff after 2 minutes inactivity
A.OFF 5' Automatic shutoff after 5 minutes inactivity

#### 4.3.8 Peripheral unit (8th menu option / see overview and notes in Section 4.1)

Peripheral devices can only be connected if the balance has been equipped with an RS232C interface. The balance automatically saves the appropriate settings (Sections 4.3.9 – 4.3.13) for every peripheral device.



Printer Connected to a printer.

Host Connection to any desired peripheral device.

Aux. display Connection of an optional auxiliary display unit (communications

parameters cannot be selected).

#### 4.3.9 Send mode (9th menu option / see overview and notes in Section 4.1)

**Note:** This menu option is only available if the "Host" setting was selected in the 7th menu option (Peripheral unit)! It specifies how a value is transferred to a peripheral device.



S. oFF Send mode switched off.

S. Stb The next possible stable value will be transferred after the « > » key

has been pressed.

S. Cont All values are transferred automatically.
S. Auto Only stable values are transferred automatically.

S. All The current value is transferred after the « > » key has been

pressed.

#### 4.3.10 Send format (10th menu option / see overview and notes in Section 4.1)

"S. SICS":

Note: This menu option is only available if the "S. oFF" setting was not selected in the 9th menu option ("Send mode")!



It sets the data transfer format.

The MT-SICS data transfer formats are used. Please refer to the "Reference Manual MT-SICS B-S/L/L-S balances 11780447".

available from your METTLER TOLEDO dealer or downloaded from

the Internet (www.mt.com/sics-classic).

More Information please find in the Section 6.3.

"S. PM"\*: The following PM balance data transfer formats are used:


S. Stb: பபபபப1.67890பg

S. Cont: Sulul 1.67890ug SDulul 1.39110ug

S. Auto: Suuuu1.67890ug S. All: uuuuu1.67890ug

\_D\_\_\_1.39110\_a

#### 4.3.11 Baud rate (11th menu option / see overview and notes in Section 4.1)

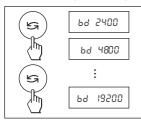

**Note:** This menu option is only available if the "Printer" or "Host" setting was selected in the 8th menu option (Peripheral unit)!


The baud rate (data transfer rate) determines the speed of transmission via the serial interface. The unit is the baud (bd) = 1 bit/second.

The following settings are available: 600 bd, 1200 bd, 2400 bd, 4800 bd, 9600 bd and 19200 bd.

For problem-free data transmission the sending and receiving devices must be set at the same value.

#### 4.3.12 Bit/Parity (12th menu option / see overview and notes in Section 4.1)

**Note:** This menu option is only available if the "Printer" or "Host" setting was selected in the 8th menu option (Peripheral unit)! It sets the character format for the peripheral device connected to the balance.

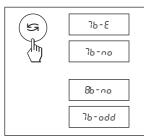

7b-E 7 data bits/even parity
7b-no 7 data bits/no parity
8b-no 8 data bits/no parity
7b-odd 7 data bits/odd parity

#### 4.3.13 Handshake (13th menu option / see overview and notes in Section 4.1)

**Note:** This menu option is only available if the "Printer" or "Host" setting was selected in the 8th menu option (Peripheral unit)! This function is used to select the data transfer mode to suit different serial devices.



HS oFF No handshake


HS SoFt Software handshake (XON/XOFF)
HS HArd Hardware handshake (RTS/CTS)

<sup>\*</sup> unidirectional, no MT-SICS commands are accepted.

#### 5 Functions

Settings and values saved under a given function are retained until they are replaced or another function is selected. The «C» key can be used to cancel the procedure currently in progress.

#### 5.1 Recall weight / Recall weight value

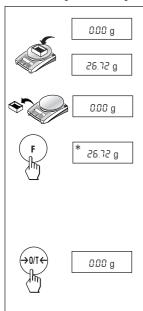

#### Requirement

The function "rEcALL" must be activated in the menu (Section 4).

- Put weight on balance.
   Display shows weight value and stores stable value.
- → Remove weight.
  When the weight is removed the Display shows zero.
- → Press the «F» key briefly

The display **shows** last stored stable weight value **for 5 seconds** together with asterisk (\*) symbol. After 5 seconds or by pressing the «**F**» key briefly, the display goes back to zero. This can be repeated unlimited times. Every recalled value is marked with the asterisk (\*) icon.


#### Delete last weight value

As soon a new stable weight value is displayed, the old value becomes replaced by the new weight value.

→ When pressing the «→0/T←» key briefly, the stored value is set to 0 and normal tare is executed.

Note: If the power is switched off, the stored value is permanently lost.

#### 5.2 Piece counting

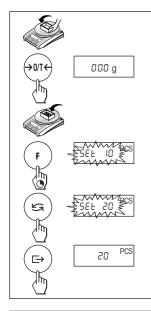


#### Requirement

The function "F count" must be activated in the menu (Section 4).

→ Place empty container on the balance and tare by briefly pressing the «→0/T←» key.

Setting the reference: a reference weight must first be entered for piece counting:


- → Add a number of reference pieces to container. Possible numbers are 5, 10, 20, 50, 100 and "no" (piece counting deactivates).
  - Note that the minimum weight = 10d (d: display increment), and the minimum unit weight = 1d!
- → Hold the «F» key down until "SEt ... PCS" is displayed.
- → Repeatedly press the «S» key until the display equals the number of reference pieces entered.
- → Confirm the number of reference pieces with the «□→» key or automatic acceptance after 7 seconds. The current number of pieces (PCS = pieces) is displayed.

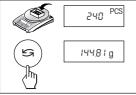

#### Switching between piece count and weight display

- → Place the items to be counted in the container. The number of pieces is displayed.
- → Press the «与» key. The weight is displayed (in unit 1, and if the key is pressed again, in unit 2, provided this function is activated).
- → Return to the piece count display by pressing the «S» key again.

#### 5.3 Percent weighing

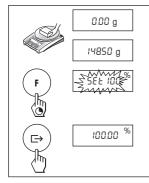


#### Requirement

The function "F 100 %" must be activated in the menu (Section 4).

#### Set target weight

- → Target weight (Reference weight, which corresponds to 100 %) in centre of pan.
  - Note that the minimum weight = 10d (d: display increment).
- → Hold the «F» key down until "SEt 100 %" is displayed.
- → Press the «与» key to select "SEt 100 %" or "SEt no %" (Percent weighing deactivated).
- → The «□→» key can be used briefly to confirm or automatic acceptance after 7 seconds.

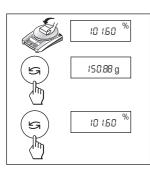

#### Switching between percent weighing and weight display

- → Place weighing sample in centre of pan.
  The weight of the sample is displayed as a percentage of the target weight.
- → Press the «S» kev. The weight is displayed.
- → Return to display in percent: pressing the «与» key again.

#### 5.4 Switching weight units



#### Requirement

Different weight units must be activated in the menu for unit 1 and unit 2 (Section 4).

→ The «≤» key can be used at any time to toggle between the two weighing units selected in the menu ("UNIT 1" and "UNIT 2").

#### Note

 Switching between weight units may be blocked with certified balances, depending on national weights and measures legislation.

#### 6 Technical data, options, optional equipment

#### 6.1 Technical data

#### Standard equipment of PL-L balances

AC adapter to national codes.
 100–240 VAC/50–60 Hz, 0.3 A
 12 VDC, 0.84 A

Balance power input 6-14,5VAC, 50/60Hz, 4VA or 7-20VDC, 4W

· All models can weigh below balance

#### **Materials**

• Housing: plastic (ABS/PC)

• Weighing pan: 18/10 chromium-nickel steel

#### **Batteries**

 4 x AA (LR6) 1.5 V alkali-manganese, typical 20 h (with 2.9 Ah capacity)

#### **Protection**

- · Protected against dust and water
- · Pollution degree: 2
- Installation category: class II
- EMC: see declaration of conformity

#### **Ambient conditions**

The technical data are valid under the following ambient conditions:

Ambient temperature 10 °C ... 30 °C

Relative humidity 10 % to 80 % at 31 °C,

linear decreasing to 50 % at 40 °C noncondensing


Operability is assured at ambient temperatures between 5 and 40  $^{\circ}\text{C}.$ 

	PL202-L	PL402-L	PL602-L
Max. capacity	210 g	410 g	610 g
Readability	0.01 g	0.01 g	0.01 g
Repeatability (sd)	0.01 g	0.01 g	0.01 g
Linearity	0.02 g	0.03 g	0.03 g
Sensitivity temperature drift (10 °C 30 °C)	±10 ppm/°C	±10 ppm/°C	±10 ppm/°C
Settling time, typical	3 s	3 s	3 s
Adjustment weight external (optional)	200 g	200 g	500 g
Interface RS232C	optional	optional	optional
External dimensions of balance (W/D/H)	194x225x67 mm	194x225x67 mm	194x225x67 mm
External dimensions of packaging (W/D/H)	350x275x140 mm	350x275x140 mm	350x275x140 mm
Weighing pan	ø 120 mm	ø 120 mm	ø 120 mm
Net weight (with packaging)	1.0 (2.1) kg	1.0 (2.1) kg	1.0 (2.1) kg
Level indicator	yes	yes	yes
Number of leveling screws	2	2	2
Certified balance available	no	no	yes

	PL601-L	PL601-L	PL1001-L	PL2001-L
Max. capacity	610 g	610 g	1100 g	2100 g
Readability	0.1 g	0.1 g	0.1 g	0.1 g
Repeatability (sd)	0.1 g	0.1 g	0.1 g	0.1 g
Linearity	0.2 g	0.2 g	0.2 g	0.2 g
Sensitivity temperature driff (10 °C 30 °C)	±10 ppm/°C	±10 ppm/°C	±10 ppm/°C	±10 ppm/°C
Settling time, typical	3 s	3 s	3 s	3 s
Adjustment weight external (optional)	500 g	500 g	1000 g	2000 g
Interface RS232C	optional	optional	optional	optional
External dimensions of balance (W/D/H) in mm	194x225x67	194x225x67	194x225x67	194x225x67
External dimensions of packaging (W/D/H) in mm	350x275x140	350x275x140	350x275x140	350x275x140
Weighing pan	ø 160 mm	ø 160 mm	ø 160 mm	ø 160 mm
Net weight (with packaging)	1.3 (2.3) kg	1.3 (2.3) kg	1.3 (2.3) kg	1.3 (2.3) kg
Level indicator	no	yes	no	yes
Number of leveling screws	_	4	_	4
Certified balance available	no	yes	no	no

	PL4001-L	PL6001-L	PL6000-L	PL6000-L
Max. capacity	4100 g	6100 g	6100 g	6100 g
Readability	0.1 g	0.1 g	1 g	1 g
Repeatability (sd)	0.1 g	0.1 g	1 g	1 g
Linearity	0.2 g	0.2 g	2 g	2 g
Sensitivity temperature driff (10 °C 30 °C)	±10 ppm/°C	±10 ppm/°C	±10 ppm/°C	±10 ppm/°C
Settling time, typical	3 s	3 s	2 s	2 s
Adjustment weight external (optional)	2000 g	5000 g	5000 g	5000 g
Interface RS232C	optional	optional	optional	optional
External dimensions of balance (W/D/H) in mm	194x225x67	194x225x67	194x225x67	194x225x67
External dimensions of packaging (W/D/H) in mm	350x275x140	350x275x140	350x275x140	350x275x140
Weighing pan	ø 160 mm	ø 160 mm	ø 160 mm	ø 160 mm
Net weight (with packaging)	1.3 (2.3) kg	1.3 (2.3) kg	1.3 (2.3) kg	1.3 (2.3) kg
Level indicator	yes	yes	no	yes
Number of leveling screws	4	4	_	4
Certified balance available	no	yes	no	yes

#### 6.2 RS232C interface

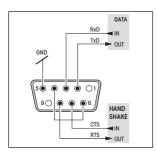

Every balance can be equipped with an optional RS232C interface for connection to a peripheral device (e.g. printer, auxiliary display or PC with a 9-pin male connector, see Section 6.4). The balance must then configured to suit the peripheral device in a menu dialog (Sections 4.3.8 - 4.3.13).

A detailed description of the available interface commands is given in the "Reference Manual MT-SICS B-S/L/L-S balances 11780447". This can be downloaded from the Internet (www.mt.com/sics-classic) and is only available in English.

The wide range of features of the PL-L balances regarding documentation of the results can be utilized by connecting to a printer, e.g. the RS-P26 or LC-P45 from METTLER TOLEDO. Printed results then make a decisive contribution to simplifying GLP/GMP-compliant work.

#### 6.3 MT-SICS Interface commands and functions

Many of the balances used have to be capable of integration in a complex computer or data acquisition system.

To enable you to integrate balances in your system in a simple manner and utilize their capabilities to the full, most balance functions are also available as appropriate commands via the data interface.

All new METTLER TOLEDO balances launched on the market support the standardized command set "METTLER TOLEDO Standard Interface Command Set" (MT-SICS). The commands available depend on the functionality of the balance.

#### Basic information on data interchange with the balance

The balance receives commands from the system and acknowledges the command with an appropriate response.

#### Command formats

Commands sent to the balance comprise one or more characters of the ASCII character set. Here, the following must be noted:

- Enter commands only in uppercase.
- The possible parameters of the command must be separated from one another and from the command name by a space (ASCII 32 dec., in this description represented as 

  ).
- The possible input for "text" is a sequence of characters of the 8-bit ASCII character set from 32 dec to 255 dec.
- Each command must be closed by C<sub>p</sub>L<sub>e</sub> (ASCII 13 dec., 10 dec.).

The characters  $C_{p}L_{pr}$ , which can be inputted using the Enter or Return key of most entry keypads, are not listed in this description, but it is essential they be included for communication with the balance.

#### Example

#### S - Send stable weight value

Command s Send the current stable net weight value.

Response SuSuWeightValueuUnit

Current stable weight value in unit actually set under unit 1.

sul Command not executable (balance is currently executing another command, e.g.

taring, or timeout as stability was not reached).

su+ Balance in overload range. su- Balance in underload range.

Example

Command s Send a stable weight value.

Response SUSUUUUU100.00Ug

The current, stable weight value is 100.00 g.

The MT-SICS commands listed below is a selected list of available commands. For additional commands and further information please refer to the Reference Manual "MT-SICS for B-S/L/L-S balances 11780447" downloadable from the Internet under www.mt.com/sics-classic.

S – Send stable weight value

Command s Send the current stable net weight value.

SI - Send value immediately

Command SI Send the current net weight value, irrespective of balance stability.

SIR - Send weight value immediately and repeat

Command SIR Send the net weight values repeatedly, irrespective of balance stability.

Z - Zero

Command **z** Zero the balance.

@ - Reset

Command @ Resets the balance to the condition found after switching on, but without a zero setting

being performed.

SR - Send weight value on weight change (Send and Repeat)

Command SR Send the current stable weight value and then send continuously the stable weight

value after every weight change.

The weight change must be at least 12.5 % of the last stable weight value,

minimum = 30d.

ST – Send stable weight after pressing □→ (transfer) key

Command ST Inquiry of actual status of the ST function.

SU - Send stable weight value with currently displayed unit

Command sv As the "s" command, but with the currently displayed unit.

#### 6.4 Optional equipment


AC adapter AC adapter universal (EU, USA, AU, UK) 100–240 VAC/50–60 Hz, 0.3 A	11120270	In-use cover All models	12102980
12 VDC, 0.84 A  AccuPac B-S Rechargeable external power source for 18 hours weighing operation independent		Printer, Application printer (LC-P45) <sup>1)</sup> Plain-paper printer, 24 characters, with additional functions (time, date, statistic, multiplier etc.)	00229119
of AC power supply	21254691	<b>Printer, Report printer (RS-P26)</b> <sup>1)</sup> Plain-paper printer, 24 characters, with	12120788
Adjustment weights  Available as OIML weights (E1, E2, F1, with calibration certificate); for further details see	11795461	additional functions (date and time).  Software 1)	
METTLER TOLEDO Weights brochure or see www.mt.com/weights		LabX direct balance (software for easy data transfer to PC)	11120340
Antitheft device Cable with lock (for all models)	00590101	Transport case For all models (without draftshield); accomodates balance,	12102982
<b>Auxiliary display <sup>1)</sup></b> Auxiliary display with switchbox	12120057	AC adapter, batteries and weights	
<b>Draft shield</b> Glass cylinder (see also weighing pan)	12102988	Weighing pan Only for models with Ø 160 mm weighing pan: Ø 120 mm weighing pan (+ pan holder + draffshield element for operation without a draff shield): necessary for use together with draft shield	12102987
Interface RS232C The interface must be fitted in the factory. Retrofi possible if carried out by a METTLER TOLEDO se	tting is only	(12102988)	

Interface cable 1)

<ul> <li>RS9–RS25: (m/f), length 2 m</li> </ul>	11101052
<ul> <li>RS9–RS9: (m/f), length 1 m</li> </ul>	11101051
<ul> <li>RS9–RS9: (m/m), length 1 m</li> </ul>	21250066
<ul> <li>RS232–USB converter cable</li> </ul>	11103691

<sup>1)</sup> RS232 interface necessary

#### 6.5 Dimensional drawing (in mm)

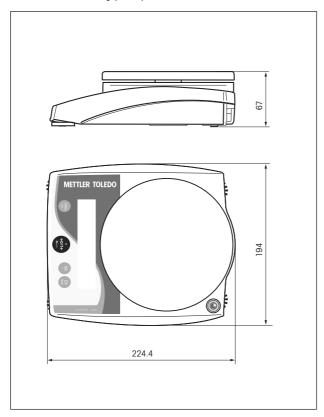

#### 7 Appendix

#### 7.1 Typical printouts from METTLER TOLEDO RS-P26 and LC-P45 printers

#### Function: Adjusting

	-
-BALANCE CALI	
METTLER TOLEDO	)
Type:	PL602-L
SNR:	1120053108
SW:	
1.0	
Weight ID:	
Weight: 200	1.00 g
External Cal.	done
Signature:	
END	

### Function: Percent weighing

I		οlρ	-	WEIGHING	G
ı	Ref.			10.008	g
ı				100.00	de
ı					
ı				60.01	g
ı				599.59	de

## Function: **Piece counting**Printout with reference weight

	PIECE	COUNTIL	٧G	
APW:		0.99	g	
Out o	f:	10		PCS
		27.00	g	
		27	PC	CS

# Function: **List**Printout of the current balance settings --- LIST OF SETTINGS ---

04.07.2006	09:55:12
METTLER TOLED Type: SNR: SW: TDNR: 7.17	PL602-L 1120053108 1.0 7.1.286.108
Weighing Para Weighing Mod Unit 1 A.Zero	
System Parame Auto off	ters:
Peripheral De P.Device Baud Bit/Parity Handshake	vices: Printer 2400 7b-even Off
P.Device Sendmode Baud Bit/Parity Handshake Soft	Host Off 9600 8b-no
END	

#### Function: Multiplier

Only possible with LC-P45. Function is triggered via the printer.

04.07.200	6 08:23:22
ID	242
SNR:	1118015657
Factor	1.65
	588.43 g
*	970.9095

## Function: Verification of the calibration (adjustment) with external weight.

Only possible with LC-P45. Function is triggered via the printer.

BALANCE TEST 04.07.2006 09:52:12
METTLER TOLEDO Type: PL602-L SNR: 1120053108 SW: 1.0
Weight ID:
Target:
External test done
Signature:
END

#### Function: Statistics

Only possible with LC-P45. Function is triggered via the printer.

04.07.200	6 10:44:07
ID	666
SNR:	1118015657
1	1100.15 g
2	1600.10 g
3	1699.95 g
n	3
х	1466.733 g
s	321.372 g
srel	21.91 %
min.	1100.15 g
max.	1699.95 g
dif.	599.80 g
	END

#### Notes

The operating instructions for the LC-P45 include a description of the functions that are triggered via that printer.

The RS-P26 prints all reports in English. This applies also to the LC-P45 reports that originate in the balance. In the case of reports triggered by the LC-P45, the following languages may be selected: German, English, French, Spanish or Italian.

#### 7.2 What if ...?

Error/Error message Cause		Rectification		
۲	Overload	→ Remove sample from weighing pan, zero again (tare).		
LJ	Underload	Check whether weighing pan is positioned properly.		
Error 1	No stability  in taring or adjusting (calibration)  when reference weight for piece counting is placed on pan	→ Wait for stability before pressing key. → Ensure more stable ambient conditions. → Remove weighing pan and clean if necessary		
Error 2	Wrong adjustment weight on pan or none at all	Place required adjustment weight in centre of pan.		
Error 3	Reference weight (Piece counting, Percent weighing) too small	→ Increase reference weight.		
Error 4	Internal fault	Contact METTLER TOLEDO customer service.		
₹0.000 g	Wrong weighing pan or pan missing or not empty	Place correct pan or empty pan on balance.		
Rbort	Adjustment aborted with the «C» key			
	No display	→ Check AC power supply.     Plug AC adapter into power supply.     → Replace batteries		

#### 7.3 Maintenance and cleaning



#### Service

Regular servicing of your balance by a service technician prolongs its working life. Ask your METTLER TOLEDO dealer for details of servicing options.

#### Cleanina

Every now and then, clean the weighing pan, draftshield element, draftshield (depending on the model) and housing of your balance using a damp cloth. Your balance is made of high-quality, durable materials and can therefore be cleaned with a standard, mild cleaning agent.



#### Please observe the following notes

- On no account use cleaning agents, which contain solvents or abrasive ingredients, as this can result in damage to the terminal overlay.
- After working with chemicals, it is advisable to wash or clean the weighing pan and the bottom
  plate (if draft shield fitted).
- Although all materials are of high quality, corrosion may occur if corrosive substances are deposited on chrome steel for an extended period of time (and if air is excluded, for example by a coating of grease).
- Ensure that no liquid comes into contact with the balance or the AC adapter!
- Never open the balance or AC adapter they contain no components, which can be cleaned, repaired or replaced by the user.
- Soiled protective covers can be replaced on all balance types (see Optional equipment).



#### Disposal

In conformance with the European Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements. Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment.

If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

Thank you for your contribution to environmental protection.

#### 7.4 Declaration of conformity

The undersigned declare on behalf of

Mettler-Toledo AG Im Lanaacher

CH-8606 Greifensee

that the balances METTLER TOLEDO PL... -L to which this declaration relates (serial number specified on the product) are in compliance with the below mentioned EEC Directives (including all amendments)

73/23/EEC Low Voltage Directive

89/336/EEC Electromagnetic compatibility and that following standards have been applied

IEC/EN61010-1:2001.

EN61326:1997+ A1:98+A2:01+A3:03 (class B)

EN61326:1997+ A1:98+A2:01+A3:03 (Minimal requirements)

for Canada, USA and Australia

CAN/CSA-C22.2 No.61010-1-04, UL Std. No.61010A-1, FCC, Part 15, class A,

AS/NZS CISPR 22, AS/NZS 61000.4.3

Balances in certified version additionally compliy with 90/384/EEC Non-automatic weighing instruments and standard EN45501.

Greifensee, 03.10.2007

Mettler-Toledo AG

Laboratory & Weighing Technologies

lengsahz

René Lenggenhager General Manager

Marcel Strotz

Manager SBU LAB Basic Weighing

To protect your METTLER TOLEDO product's future:
METTLER TOLEDO Service assures the quality, measuring accuracy
and preservation of value of all METTLER TOLEDO products for years to
come.

Please send for full details about our attractive terms of service. Thank you.



Subject to technical changes and to the availability of the accessories supplied with the instruments.