Operating Instructions

Excellence Plus Precision Balances

XP Models - Part 1

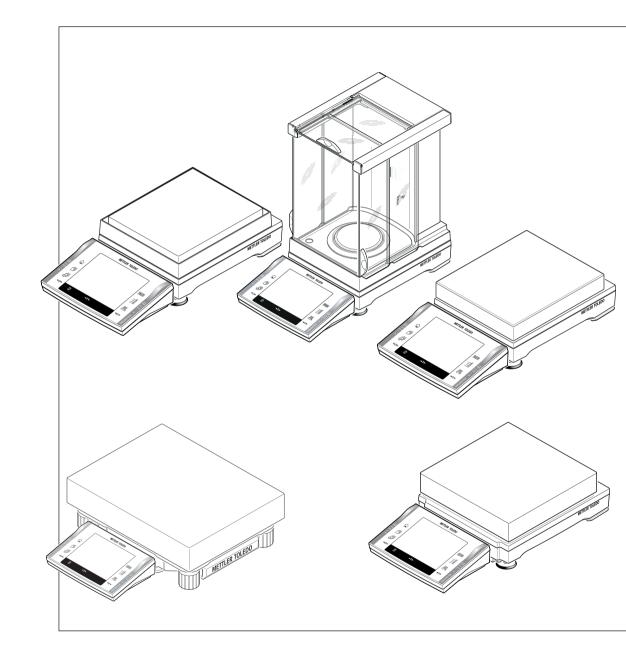




Table of contents

1	Introduction	n		5
		1.1	Conventions and Symbols Used in These Operating Instructions	6
2	Safety Info	mation		7
		2.1	Definition of Signal Warnings and Symbols	7
		2.2	Product Specific Safety Notes	7
3	Overview X	P Precisio	n Balances	9
		3.1	Overview "S" and "M" Weighing Platform	9
		3.2	Overview "L" Weighing Platform	11
4	Setting up t	the Balanc	e	13
		4.1 4.1.1	Unpacking and Checking the Delivered Items Unpacking the Draft Shield "S" Weighing Platform (Model-dependent)	13 13
		4.2	Scope of Delivery	13
		4.3	Selecting a Location	15
		4.4	Assembling the Balance	15
		4.4.1	Installing the Balance Cable in the Terminal "S" and "M" Weighing Plat- form	15
		4.4.2 4.4.3	Placing the Terminal on the "S" and "M" Weighing Platform Fastening the Terminal to the Terminal Support with Screws "S" and "M" Weighing Platform	16 17
		4.4.4	Placing the Terminal on the "L" Weighing Platform	17
		4.4.5	Installing the Draft Shield and Weighing Pan "S" and "M" Weighing Plat- form	19
		4.5	Power Supply	20
		4.5.1 4.5.2	Power Supply for the "S" and "M" Weighing Platform Power Supply for the "L" Weighing Platform	20 21
		4.6	Optimise the Readability of the Terminal and Location of the Terminal	22
		4.6.1	Setting the Reading Angle	22
		4.6.2	Removing the Terminal and Placing it Within Range of the Weighing Plat- form	22
		4.6.2.1 4.6.2.2	Setting up the Terminal Separately "S" and "M" Weighing Platform Setting up the Terminal Separately "L" Weighing Platform	22 23
		4.7 4.7.1 4.7.2	Transporting the Balance Transporting Over Short Distances "S" and "M" Weighing Platform Transporting Over Short Distances "L" Weighing Platform	23 23 24
		4.7.3 4.8	Transporting Over Long Distances	24
		4.8 4.8.1 4.8.2	Below-the-Balance Weighing Below-the-Balance Weighing "S" and "M" Weighing Platform Below-the-Balance Weighing "L" Weighing Platform	25 25 25
5	First Steps			27
•		5.1	Switching On / Off	27
		5.2	Leveling the Balance	27
		5.2.1	Leveling the Balance 10 mg, 0.1 g and 1 g "S" and "M" Weighing Plat- form	28
		5.2.2 5.2.3	Leveling the Balance 1 mg "S" and "M" Weighing Platform Leveling the Balance "L" Weighing Platform	29 29

6	Maintenan	ce		31
		6.1	Cleaning	31
		6.2	Cleaning the Draft Shield (0.1 mg and 1 mg Models)	31
		6.3	Disposal	32
7	Technical I	Data		33
		7.1	General Data "S" and "M" Weighing Platform	33
		7.2	Explanatory Notes for the METTLER TOLEDO AC Adapter	33
		7.3	General Data "L" Weighing Platform	34
		7.4	Model-specif Data	35
		7.4.1 7.4.2 7.4.3 7.4.4 7.4.5 7.4.6	Balances with Readability of 0.1 mg, "S" Platform with Draft Shield Balances with Readability of 1 mg, "S" Platform with Draft Shield Balances with Readability of 10 mg, "S" Platform with Draft Shield Element Balances with Readability of 0.1 g, "S" Platform Balances with Readability of 10 mg / 0.1 g / 1 g, "M" Platform Balances with Readability of 0.1 g / 1 g, "L" Platform	35 36 38 41 43 47
		7.5 7.5.1 7.5.2 7.5.3 7.5.4 7.5.5 7.5.6	Dimensions Balances with Readability of 0.1 mg, "S" Platform with Draft Shield Balances with Readability of 1 mg, "S" Platform with Draft Shield Balances with Readability of 10 mg, "S" Platform with Draft Shield Element Balances with Readability of 0.1 g, "S" Platform Balances with Readability of 10 mg / 0.1 g / 1 g, "M" Platform Balances with Readability of 0.1 g / 1 g, "L" Platform	50 50 51 52 53 54 56
		7.6 7.6.1 7.6.2	Interfaces Specifications of RS232C Specifications of "Aux" Connection	57 57 57
8	Accessorie	s and Sp	are Parts	58
		8.1	Accessories	58
		8.2	Spare Parts	68
9	Appendix			70
		9.1	MT-SICS Interface Commands and Functions	70
		9.2	Procedure for Certified Balances	70
10	Index			72

1 Introduction

Thank you for choosing a METTLER TOLEDO balance.

The balances of the XP line combine a large number of weighing and adjustment possibilities with exceptionally convenient operation.

In this chapter you will be given basic information about your balance. Please read right through this chapter carefully even if you already have experience with METTLER TOLEDO balances. Please pay special attention to the safety warnings!

The different models have different characteristics regarding equipment and performance. Special notes in the text indicate where this makes a difference to operation.

The XP line comprises a range of balances which differ from each other in relation to their weighing range and resolution.

The following features are common to all models of the XP lines:

- Fully automatic adjustment "ProFACT" using internal weights.
- Built-in level sensor, illuminated level indicator and Leveling Assistant for fast and easy leveling.
- Built-in applications for normal weighing, statistics, formulation, piece counting, percent weighing, density, dynamic weighing, differential weighing and LabX Client.
- Integral RS232C interface.
- Slot for second interface (optional).
- Touch-sensitive graphics terminal ("Touch Screen") with color display.
- Two programmable sensors for hands-off operation ("SmartSens") to speed up frequently recurring tasks.

A brief word about standards, guidelines, and methods of quality assurance: The balances comply with usual standards and guidelines. They support standard procedures, specifications, working methods, and reports according to **GLP** (**G**ood Laboratory **P**ractice). In this connection, records of working procedures and adjustments become very important; for this purpose we recommend you to use a printer from the METTLER TOLEDO range, since these are optimally adapted to your balance. The balances conform to the applicable standards and guidelines and possess a EC declaration of conformity. METTLER TOLEDO is certified as manufacturer according to ISO 9001 and ISO 14001.

The Operating Instructions for the XP balances consist of 3 separate documents, whose contents are listed bellow.

Part 1, This Document Contents

- Introduction
- Safety Information
- Setting up the Balance
- Leveling the Balance
- Cleaning and Service
- Technical Data
- Interface commands and MT-SICS functions
- Accessories
- Spare Parts

Part 2, Separate Document

Contents: Terminal, System and Applications

- Basic Principles for Using the Terminal and the Firmware
- System Settings
- User-specific Settings
- Applications
- Firmware (Software) Updates
- Error and Status Messages
- Conversion Table for Weight Units
- Recommended Printer Settings

Part 3, Separate Document Contents: Adjustments and Tests

- Adjustments
- Tests

Finding More Information

Internet http://www.mt.com/excellence

1.1 Conventions and Symbols Used in These Operating Instructions

The following conventions apply to the operating instructions: Part 1, Part 2 and Part 3.

Key designations are indicated by a picture or text in double angular parentheses (e.g. «==» or «On/Off»).



This symbol indicates press key briefly (less than 1.5 s).



This symbol indicates press and hold key down (longer than 1.5 s).

These symbols indicate an instruction:

- prerequisites
- 1 steps
- 2 ...
- → results

2 Safety Information


2.1 Definition of Signal Warnings and Symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the balance, malfunctions and false results.

Signal Words

WARNING	for a hazardous situation with medium risk, possibly resulting in severe injuries or death if not avoided.
CAUTION	for a hazardous situation with low risk, resulting in damaged to the device or the property or in losing of data or minor or medium injuries if not avoid- ed.
Attention	(no symbol) for important information about the product.
Note	(no symbol) for useful information about the product.

Warning Symbols

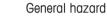



Electrical shock

2.2 Product Specific Safety Notes

Always operate and use your balance only in accordance with the Operating Instructions Part 1, Part 2 and Part 3.

The instructions for setting up your new balance must be strictly observed.

If the instrument is not used according to the manufacturer's Operating Instructions (Part 1, Part 2 and Part 3), protection of the instrument may be impaired.

Intended Use

Your balance is used for weighing. Use the balance exclusively for this purpose. Any other type of use and operation beyond the limits of technical specifications without written consent from Mettler-Toledo AG, is considered as not intended.



It is not permitted to use the instrument in explosive atmosphere of gases, steam, fog, dust and flammable dust (hazardous environments).

CAUTION



Damage of Device

- For use only in dry interior rooms.

 Do not use sharply pointed objects to operate the keyboard! Although your balance is very ruggedly constructed, it is nevertheless a precision instrument. Treat it with corresponding care.

- Do not open the balance: It does not contain any parts which can be maintained, repaired, or replaced by the user. If you ever have problems with your balance, contact your METTLER TOLEDO dealer.

 Use only balance accessories and peripheral devices from METTLER TOLEDO; they are optimally adapted to your balance.

CAUTION




Damage of Device

Use only the original universal AC adapter delivered with your balance, and check that the voltage printed on it is the same as your local power supply voltage. Only plug the adapter into a socket which is grounded.


3 Overview XP Precision Balances

3.1 Overview "S" and "M" Weighing Platform

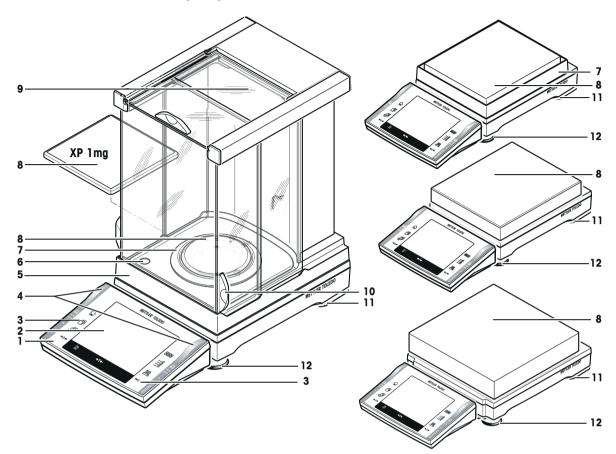

Front View S and M Platform


1	Terminal (details see Operating Instruction – Part 2)	2	Display (Touch-sensitive "Touch Screen")
3	Operating keys	4	SmartSens sensors
5	Type name	6	Level indicator / level sensor
7	Draft shield element	8	Weighing pan
9	Glass draft shield	10	Handle for operation of the draft-shield door
11	Safety feet (10 mg, 0.1 g + 1 g models)	12	Foot screws

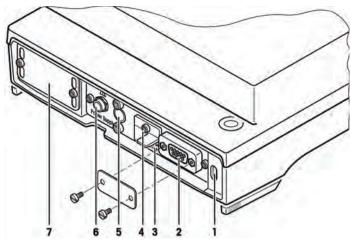

Rear View S and M Platform


1	Fastening point for anti-theft device	2	RS232C serial interface
3	Aux 1 (connection for "ErgoSens", hand- or foot-switch)		Aux 2 (connection for "ErgoSens", hand- or foot-switch)
5	Fastening for auxiliary display stand or termi- nal stand (optional)	6	Socket for AC adapter
7	Slot for second interface (optional)		

3.2 Overview "L" Weighing Platform

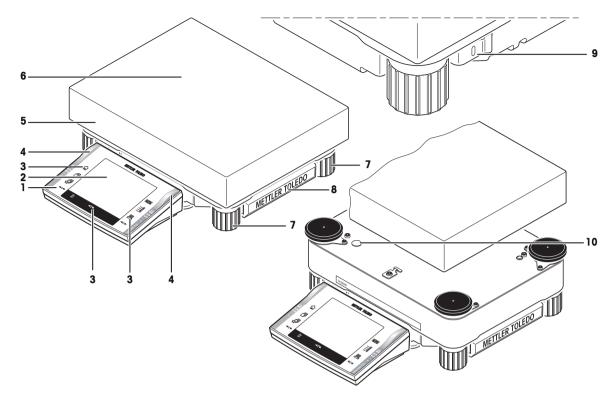

Top View L Platform

1	Terminal (details see Operating Instruction – Part 2)	2	Display (Touch-sensitive "Touch Screen")
3	Operating keys	4	SmartSens sensors
5	Type name	6	Weighing pan
7	Foot screws	8	Cover
9	Fastening point for anti-theft device	10	Level indicator / level sensor

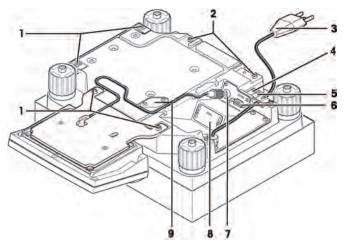

Bottom View L Platform

1	Points of attachment for terminal or cover	2	Fastening for terminal stand (optional)
3	Power cable	4	Aux 1 (connection for "ErgoSens", hand- or foot-switch)
5	Aux 2 (connection for "ErgoSens", hand- or foot-switch)	6	RS232C serial interface
7	Connector for terminal cable	8	Slot for second interface (optional)
9	Cover plate for below-the-balance weighing (hook optional)		

4 Setting up the Balance

This chapter describes how to unpack your new balance, set it up and prepare it for operation. On completion of the steps described in this chapter, your balance is ready for operation.



CAUTION

Electrical shock

The balance must be disconnected from the power supply when carrying out all setup and mounting work.

4.1 Unpacking and Checking the Delivered Items


Open the packaging and carefully remove all components.


4.1.1 Unpacking the Draft Shield "S" Weighing Platform (Model-dependent)

- 1 Place the draft shield on a clean surface.
- 2 Turn so the cover (1) is vertically on top.
- 3 Lift the carton (2) off over the handle and the pull the carton off toward the back.
- 4 Hold the glass panels firmly so they do not get pulled off and fall to the floor.
- 5 Close the cover (1) again.
- 6 Push all the glass panels as far as they will go to the back.

Balance with 0.1 mg readability

- 1 Place the ring seal (3) into the draft shield base (4) from above.
- 2 Push the complete ring seal (3) through the opening in the draft shield base.
- 3 Pull the top edge upward through the opening until the edge lies flat and tight all the way round.
- 4 Check by running a finger all the way round that the seal ring (3) is fitted tightly into the opening in the draft shield base (4).
- 5 Insert the bottom plate (5).

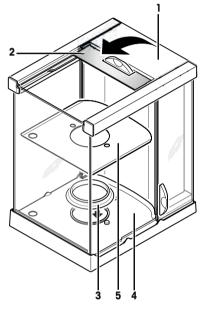

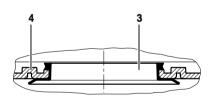

Balance with 1 mg readability

Insert the bottom plate (5).

4.2 Scope of Delivery

The standard scope of delivery contains the following items:

Components		S platform	1		
		0.1 mg	1 mg	10 mg	0.1 g
Weighing plafform		 ✓ 	\checkmark	1	1
	RS232C interface	 ✓ 	1	1	1
	Slot for second interface (optional)	1	\checkmark	1	1

Components			S platform	1		
	Feedthroughs for below-the-bala weighing	nce	<i>✓</i>	1	√	1
	Feedthroughs for antitheft device		 ✓ 	1	 ✓ 	\checkmark
AC adapter			1	1	 ✓ 	1
Country-specific pow	ver cable		1	1		1
Terminal support			1	1		1
Terminal with protect	tive cover		 ✓ 	 ✓ 	 ✓ 	1
Draft shield with bott	om plate		 ✓ 	1	_	_
Draft shield element				_		_
Ring seal				_		_
Protective cover for v	veighing platform		-	_		 ✓
Weighing pan suppo			_	1		
Weighing pan	ø 90 mm			_	-	_
troiginig pari	127 x 127 mm		• •	 ✓ 	_	
	170 x 205 mm			•		
	190 x 223 mm		_		✓	
	237 x 237 mm		_	_		1
	280 x 360 mm		_			
Operating Instruction	s Part 1 (this document), Part 2 a	nd Part 3				
Production certificate				· /		<u> </u>
CE declaration of cor						
		M platfor	•	•	-	-
Components		10 mg	0.1 g	1 g	L platform 0.1 g	1 g
Weighing platform			v.r g	. <u> </u>	V.1 g	
0 01	RS232C interface					<u> </u>
	Slot for second interface (optional)	✓ ✓	✓ ✓	✓ ✓	✓ ✓	✓ ✓
	Feedthroughs for below-the-bal- ance weighing	1	~	\checkmark	_	-
	Ready for below-the-balance weighing (hook optional)	_	-	_	1	1
	Feedthroughs for antitheft	1	1	1	1	\checkmark
	device					
AC adapter	device	✓	✓	1	_	_
AC adapter Country-specific pow		\ \ \	\ \ \	\ \	-	-
•		-			-	-
Country-specific pow	/er cable	1	1	1	-	-
Country-specific pow Terminal support	ver cable	<i>J</i>	\ \ \	√ √	✓	\checkmark
Country-specific pow Terminal support Terminal with protect	ver cable	<i>J</i>	\ \ \	√ √	✓	\checkmark
Country-specific pow Terminal support Terminal with protect Draft shield with bott Draft shield element Ring seal	ver cable tive cover om plate	✓ ✓ ✓ −	✓ ✓ ✓ −	√ √	✓	✓ ✓ ─
Country-specific pow Terminal support Terminal with protect Draft shield with bott Draft shield element	ver cable tive cover om plate	✓ ✓ ✓ − −	✓ ✓ ✓ − −	✓ ✓ ✓ – –	✓ ✓ − −	✓ ✓ ─

Setting up the Balance	15

Components		M platform	n *		L platform	1
Weighing pan	ø 90 mm	-	-	-	-	-
	127 x 127 mm	-	-	-	-	-
	170 x 205 mm	-	-	-	-	-
	190 x 223 mm	-	-	-	-	_
	237 x 237 mm	 ✓ 	1	1	-	-
	280 x 360 mm	-	-	-	1	1
Operating Instruct and Part 3	ons Part 1 (this document), Part 2	1	1	1	1	1
Production certific	ate	 ✓ 	1	1	1	1
CE declaration of	conformity		1			

* The precision balances with M platform are no longer available.

4.3 Selecting a Location

Choose a position which is stable, free from vibration, and as nearly horizontal as possible. The supporting surface must be able to bear the weight of the fully loaded balance safely.

Avoid the following:

- Direct sunlight
- Draft (e.g. from fans or air conditioning)
- Excessive fluctuations in temperature.

Further information can be found in Weighing the Right Way.

Observe ambient conditions. See Technical Data (page 33).

Note

If the balance is not horizontal from the beginning, it will have to be leveled during initial operation. **See** Leveling the Balance (page 27).


4.4 Assembling the Balance


The terminal is identical for all XP precision balances. The size of the weighing pan depends on the readability and maximum capacity of the balance.


4.4.1 Installing the Balance Cable in the Terminal "S" and "M" Weighing Platform

- Ensure that a soft, clean surface is used so as not to damage the terminal surface.
- 1 Place the terminal on the operating surface.
- 2 Open the housing by pressing the 2 buttons (1) for adjusting the terminal and swiveling the housing base upwards.
- 3 Pull the cable with the retaining ring (2) through the opening on the housing base.
- 4 Return the terminal to its normal position.




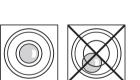


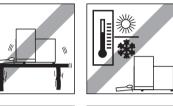

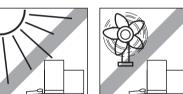

- 5 Open it so that the cable can be accessed.
- 6 Insert the cable (3) into the top housing.
- 7 Close both parts of the housing until the retaining ring (2) is positioned in the cable opening of the housing base.
- 8 Place the retaining ring (2) behind the two guideways and check that it is secure (tension relief).
- 9 Before closing the housing, check that the plug is inserted correctly into the terminal plug-in connection.
- 10 Now close the housing by pressing the two buttons (1) for adjusting the terminal until the housing base engages in the top housing.

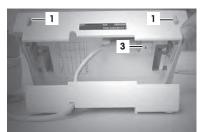

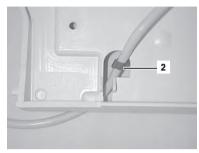




4.4.2 Placing the Terminal on the "S" and "M" Weighing Platform

 Slide the terminal support (4) under the front of the weighing platform until you feel it snap into place.
 Note

For this process, the weighing platform must not be lifted but should ideally be placed at the weighing location.

- 2 Push the terminal support in "horizontally" with it lying on the table.
- 3 Place the terminal in the center of the terminal support.
- 4 Slide it towards the weighing platform until the terminal tilts slightly downwards.

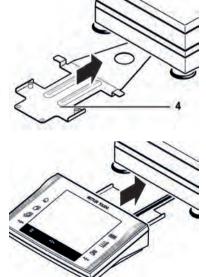

Attention

The balance and the terminal are not fastened together by the terminal support! When transporting by hand, always hold the balance and the terminal firmly, **see** Transporting the Balance (page 23).

Note

You can also place the terminal free of the terminal support anywhere around the weighing platform where the lenght of the cable allows.

4.4.3 Fastening the Terminal to the Terminal Support with Screws "S" and "M" Weighing Platform

If you change the weighing site often, we recommend fastening the terminal to the terminal support with screws.

- 1 Pull the terminal, along with the terminal support, about 2 inches (5 cm) away from the weighing platform.
- 2 Pull the terminal cable as far as possible in the direction of the terminal.
- Open the terminal by pressing the 2 buttons.


 Fasten the terminal to the terminal support using the 2 knurled screws (included in delivery).


- Before closing the terminal, check that the terminal plug is firmly in place.
- 1 The retaining ring for the cable must be correctly positioned when the terminal is closed.
- 2 Slide the terminal support with the fastened terminal under the weighing platform (do not lift the balance!) until you feel it snap into place.

4.4.4 Placing the Terminal on the "L" Weighing Platform

The terminal can be attached to the long or short side of the balance.

- 1 Mount the weighing pan.
- 2 Carefully turn the weighing platform over onto the weighing pan.

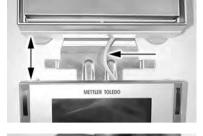

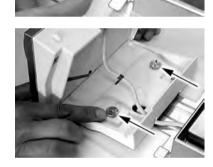






Attaching the terminal to the long side


- 1 Dismantle the cover (1) by removing the 2 screws (2).
- 2 Attach the terminal to the terminal support (5) as shown, using the screws (2) from the cover that was removed.
- 3 Insert the terminal cable (3) into the cable channel, as shown.
- 4 Screw the connector of the terminal cable into the socket (4).
- 5 Turn the balance back over into position.

2



- 1 Dismantle the cover (1) by removing the 2 screws (2).
- 2 Attach the terminal with terminal support to the points of attachment(6) with the screws (2).
- 3 Insert the terminal cable (3) into the cable channel, as shown.
- 4 Screw the connector of the terminal cable into the socket (4).
- 5 Turn the balance back over into position.

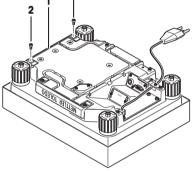

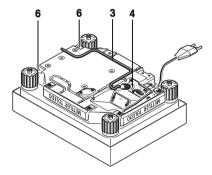

4.4.5 Installing the Draft Shield and Weighing Pan "S" and "M" Weighing Platform

Balances with readability of 0.1 mg, "S" platform with draft shield

 Place the following components on the balance in the specified order:

Attention

Push the side glass back as far as it will go. Grasp the draft shield with both hands on the bars at the top.

- Draft shield (1) with ring seal inserted, see Unpacking the Draft Shield "S" Weighing Platform (Model-dependent) (page 13).
- Bottom plate (2), if not already inserted, see Unpacking the Draft Shield "S" Weighing Platform (Model-dependent) (page 13)!
- Draft shield element (3).
- Weighing pan (4).

Note

Cleaning the draft shield, **see** Cleaning the Draft Shield (0.1 mg and 1 mg Models) (page 31).

Balances with readability of 1 mg, "S" platform with draft shield

Place the following components on the balance in the specified order:

Attention

Push the side glass back as far as it will go. Grasp the draft shield with both hands on the bars at the top.

- Draft shield (1) with ring seal inserted.
- Bottom plate (2), if not already inserted, see Unpacking the Draft Shield "S" Weighing Platform (Model-dependent) (page 13)!
- Pan support (3).
- Weighing pan (4).

Note


Cleaning the draft shield, **see** Cleaning the Draft Shield (0.1 mg and 1 mg Models) (page 31).


Balances with readability of 10 mg, "S" platform with draft shield element


- Place the following components on the balance in the specified order:
- Pan support (1).
- Weighing pan (2).
- Draft shield element (3).

Note

It is also possible to work without the draft shield element (3). However, depending on the ambient conditions, the display of the results may be slightly more unstable.

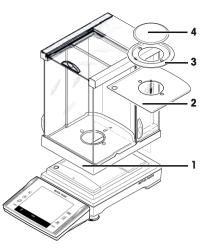


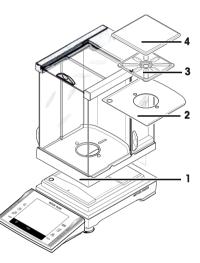

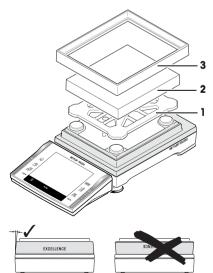

Balances with readability of 10 mg ("M" platform), 0.1 g ("S" and "M" platform) and 1 g ("M" platform)

- Place the following components on the balance in the specified order:
- Pan support (1).
- Weighing pan (2).

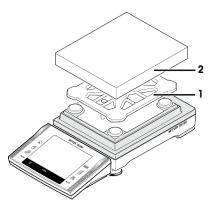

4.5 Power Supply



WARNING

Risk of electric shock


- Ensure that the AC power pack for the balance is only used in accordance with the specifications listed in chapter General Data.

- Your instrument is supplied with a 3-pin power cable with an equipment grounding conductor. Only extension cables which meet this relevant standards and also have an equipment grounding conductor may be used. Intentional disconnection of the equipment grounding conductor is prohibited.

4.5.1 Power Supply for the "S" and "M" Weighing Platform

- Your balance is delivered complete with an AC adapter and a country-specific power supply cable. The AC adapter is suitable for all power supply voltages in the range of: 100 240 VAC, 50/60 Hz, for exact specifications, see Technical Data (page 33).
- First, check whether the voltage of the power supply matches your local line voltage. If this is not the case, on no account connect the balance to the power, but contact your METTLER TOLEDO sales representative or dealer.
- Guide the cables so that they cannot become damaged and will not be in your way during your daily work! Take care that the AC adapter cannot come into contact with liquids!
- The power plug must be always accessible.
- Before operating, check all cables for damage.


- Plug the AC adapter (1) into the socket (2) in the back of your balance and into the power supply.
- After the balance has been connected to the power supply, it carries out a self test and is then ready for operation.

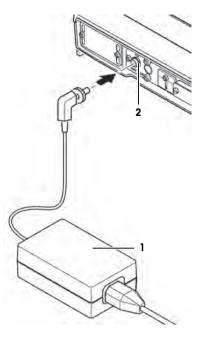

Note

If the display field remains dark, even though the power supply connection functions.

- 1 First disconnect the balance from the power supply.
- 2 Open the terminal.
- 3 Press both buttons (3) on the back of the terminal and open the upper part of the terminal.
- 4 Check that the plug for the terminal cable (4) is connected correctly inside the terminal.

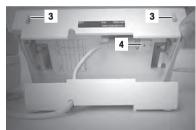

4.5.2 Power Supply for the "L" Weighing Platform

- Your balance is delivered with a country-specific power supply cable.
- First, check whether the voltage of the power supply matches your local line voltage. If this is not the case, on no account connect the balance to the power, but contact your METTLER TOLEDO sales representative or dealer.
- Guide the cables so that they cannot become damaged and will not be in your way during your daily work! Take care that the AC adapter cannot come into contact with liquids!
- The power plug must be always accessible.
- Before operating, check all cables for damage.
- Connect the balance to the power supply.
- ⇒ After the balance has been connected to the power supply, it carries out a self test and is then ready for operation.

Note

If the display field remains dark, even though the power supply connection functions.

- 1 First disconnect the balance from the power supply.
- 2 Open the terminal.
- 3 Press both buttons (1) on the back of the terminal and open the upper part of the terminal.
- 4 Check that the plug for the terminal cable (2) is connected correctly inside the terminal.

4.6 Optimise the Readability of the Terminal and Location of the Terminal

4.6.1 Setting the Reading Angle

- 1 Press in the two buttons (1) on the back of the terminal.
 - ⇒ The top of the terminal can then be pulled up or pushed down until it engages in the desired position. A total of 3 setting positions are available.
- 2 Move it in an appropriate position.

4.6.2 Removing the Terminal and Placing it Within Range of the Weighing Platform

The terminal is connected to the weighing platform with a cable. The terminal can be removed from the weighing platform and located separately to ensure the optimum setup of your workplace.


If you would like to place the terminal farther away from the balance, we recommend that you use the extension cable from our accessories range, **see** Accessories (page 58).


- 1 Switch off the balance.
- 2 Disconnect power cable.


4.6.2.1 Setting up the Terminal Separately "S" and "M" Weighing Platform

- 1 Dismantle the terminal (terminal terminal-support) by removing the 2 screws, if the terminal is mounted.
- 2 Carefully lift the terminal from the terminal support.
- 3 Pull the terminal support out of the weighing platform.
- 4 Tilt the weighing platform high on its side and carefully pull the cable out of the cable opening under the weighing platform.
- 5 Place the terminal in the desired position.

The cable can also be led out the back or the side of the weighing platform.

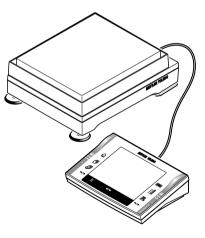

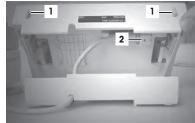


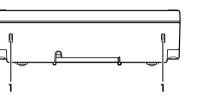

4.6.2.2 Setting up the Terminal Separately "L" Weighing Platform

- 1 Carefully turn the weighing platform over onto the weighing pan.
- 2 Carefully pull the terminal cable (3) out of the cable channel.
- 3 Remove the screws (2).
- 4 Lift the terminal with terminal support (1) off the weighing platform.
- 5 Open the housing by pressing the 2 buttons (4) for adjusting the terminal.
- 6 Unplug the cable (5) and pull it out of the housing through the opening.
- 7 Remove the 2 knurled screws (6) and detach the terminal support.
- 8 Pull the cable back through the base of the housing and plug it in.
- 9 Close the terminal and place it in the desired position.
- 10 If possible, insert the terminal cable (3) back into the cable channel.
- 11 Turn the balance back over into position.

The cable can also be led out the back or the side of the weighing platform.

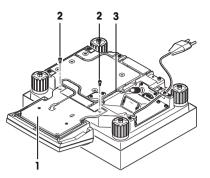





4.7 Transporting the Balance

Observe the following instructions to transport your balance to a new location.

- 1 Switch off the balance.
- 2 The balance must be disconnected from the power supply.
- 3 Remove any interface cable from the balance.

4.7.1 Transporting Over Short Distances "S" and "M" Weighing Platform

If you wish to move your balance over a short distance to a new location, proceed as follows.



CAUTION


Damage of Device

Never lift the balance by the glass draft shield, as this can cause damage! The draft shield is not fastened to the balance.

The terminal is not permanently connected to the weighing platform and therefore the balance must be carried horizontally.

 Remove the terminal from the terminal support and place the terminal on the weighing pan.

- 1 Hold the weighing platform with both hands and lift horizontally.
- 2 Carry the balance horizontally to the new location, observe the notes in chapter Selecting a Location (page 15).

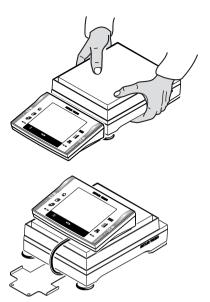

4.7.2 Transporting Over Short Distances "L" Weighing Platform




CAUTION

Damage of Device

The terminal is not permanently connected to the weighing platform and therefore the balance must be carried horizontally.

- Remove the terminal from the terminal support and place the terminal on the weighing pan.
- 1 Hold the weighing platform with both hands and lift horizontally.
- 2 Carry the balance horizontally to the new location, observe the notes in chapter Selecting a Location (page 15).

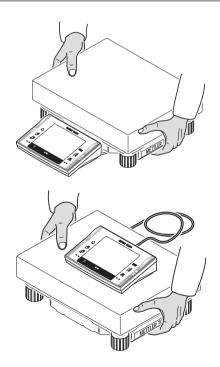

4.7.3 Transporting Over Long Distances

If you want to transport or ship your balance over long distances, or if it is not certain that the balance will be transported upright, use the complete original packaging.

4.8 Below-the-Balance Weighing

Your balance is equipped with a hanger for carrying out weighings below the work surface (weighing below the balance).

- 1 Switch off the balance.
- 2 The balance must be disconnected from the power supply.
- 3 Remove any interface cable from the balance.

4.8.1 Below-the-Balance Weighing "S" and "M" Weighing Platform

- 1 Remove the weighing pan (0.1 mg models).
- 2 Remove the draft shield element (0.1 mg and 10 mg models [S platform only]).
- 3 Remove the weighing pan and pan support.

4 Attention

Models with a glass draft shield: Carefully lift the draft shield from the weighing platform and put it aside.

- 5 Lift the terminal from the terminal support.
- 6 Remove the 2 screws between terminal and terminal support, **see** Fastening the Terminal to the Terminal Support with Screws "S" and "M" Weighing Platform (page 17).
- 7 Place the terminal to the side of the weighing platform.
- 8 Remove the terminal support.
- 9 Tilt the weighing platform backwards until the cover plate (2) is visible.

Attention

Do not place the weighing platform on the location bolt for the pan support (0.1 mg and 1 mg models).

- 10 Remove the 2 screws (1) and the cover plate (2).
 - \Rightarrow The hanger is now accessible.
- 11 Return the balance to its normal position and simply reinstall all components in the reverse order.


Note


For below-the-balance weighing with the M model (XP16001M, XP16001MDR, XP20001M and XP20000M) you will need hook 11132565 from the accessories range.

4.8.2 Below-the-Balance Weighing "L" Weighing Platform


Note

For below-the-balance weighing you will need hook 11132565 from the accessories range.

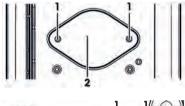

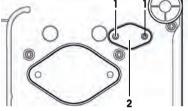

- 1 Carefully turn the weighing platform over onto the weighing pan.
- 2 Remove the 2 screws (1) and the cover plate (2).
 - \Rightarrow The hanger is now accessible.
- 3 Fasten the hooks (optional) with screws.
- 4 Return the balance to its normal position and simply reinstall all components in the reverse order.

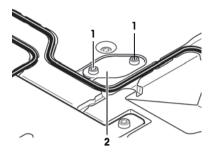


5 First Steps

5.1 Switching On / Off

Switching On

- Press «On/Off».
- \Rightarrow The display appears.

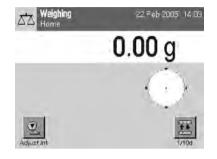

Note

If the balance has not been set up exactly horizontally, a warning text will appear shortly after the balance is turned on, prompting you to level the balance.

Switching Off

- Press «On/Off» until "Off" appears in the display.

Note

Do not disconnect the balance from the power supply except if you will not be using the balance for an extended period.

5.2 Leveling the Balance

Your balance has a built-in levelcontrol, that constantly checks the correct horizontal alignment.

When the levelcontrol detects that the level is incorrect, a warning will appear and a warning beep will sound. A status icon will also appear in the upper right-hand corner of the display.

- To start the Leveling Assistent, tap «Info».
 - ⇒ The Leveling Assistant will guide you step by step through the leveling process.






On off







5.2.1 Leveling the Balance 10 mg, 0.1 g and 1 g "S" and "M" Weighing Platform


- 1 The Leveling Assistant will prompt you to loosen the clamps for the safety feet.
- 2 Remove the clamps (1) for the safety feet by turning them outwards.
- 3 Turn the clamps (1) outwards as far as they will go (~ 90°), so that the safety feet can move freely.
- 4 Tap **«Forward >**» after you have loosened the clamps for the safety feet.

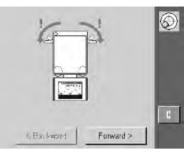

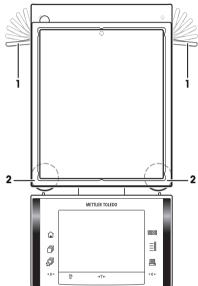

Please press the appropriate button to show the laveling direction.

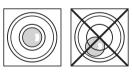

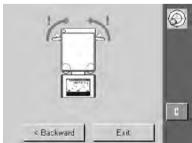

- 1 Watch the levelcontrol located of your balance and press the appropriate button of the current position.
 - ⇒ The Leveling Assistant will now use red arrows to indicate the direction in which you must turn the two footscrews (2).
- 2 Screw the footscrew until the air bubble is in the inner circle.
- 3 Tap «Forward >».
- 4 The Leveling Assistant will now prompt you to lock the clamps for the safety feet in place.
- 5 Secure the safety feet by turning the clamps inward as far as they will go.
- 6 Tap «**Exit**».
 - \Rightarrow A message appears that advises you to adjust the balance.
- 7 Confirm with «OK».
- ⇒ Status icon will no longer appear and balance returns to normal operation.

5.2.2 Leveling the Balance 1 mg "S" and "M" Weighing Platform


- 1 Watch the levelcontrol located of your balance and press the appropriate button of the current position.
 - ⇒ The Leveling Assistant will now use red arrows to indicate the direction in which you must turn the two footscrews.
- 2 Screw the footscrew until the air bubble is in the inner circle.
- 3 Tap «Exit».


 \Rightarrow A message appears that advises you to adjust the balance.


- 4 Confirm with «OK».
- ⇒ Status icon will no longer appear and balance returns to normal operation.


5.2.3 Leveling the Balance "L" Weighing Platform


- 1 Touch the button that matches your site and your configuration.
- 2 Remove the weighing pan so you can monitor the level indicator.
- 3 Tap «Forward >».
- 4 Turn the leveling foot all the way up, in the direction of the red arrow.
- 5 Tap «Forward >».
- 6 Watch the levelcontrol located of your balance and press the appropriate button of the current position.
- 7 The Leveling Assistant will now use red arrows to indicate the direction in which you must turn the footscrews.
- 8 Screw the footscrew until the air bubble is in the inner circle.
- 9 Tap «Forward >».

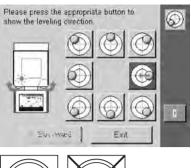


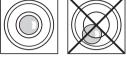

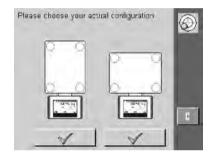


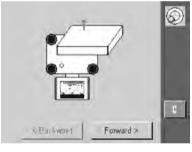


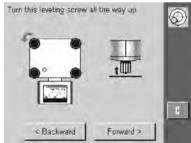


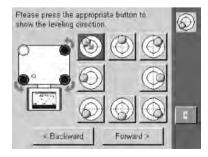

- 1 Touch the button that matches your site and your configuration.
- 2 Remove the weighing pan so you can monitor the level indicator.
- 3 Tap «Forward >».
- 4 Turn the leveling foot all the way up, in the direction of the red arrow.
- 5 Tap «Forward >».
- 6 Watch the levelcontrol located of your balance and press the appropriate button of the current position.
- 7 The Leveling Assistant will now use red arrows to indicate the direction in which you must turn the footscrews.
- 8 Screw the footscrew until the air bubble is in the inner circle.
- 9 Tap «Forward >».
- 1 Unscrew the footscrew until it **lightly** touches the supporting surface.
- 2 Tap «Forward >».
- 3 Tap «Exit».
 - \Rightarrow A message appears that advises you to adjust the balance.
- 4 Confirm with «OK».
- ⇒ Status icon will no longer appear and balance returns to normal operation.

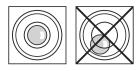

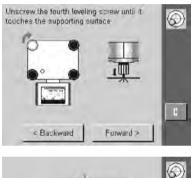

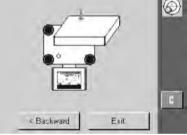

6 Maintenance

6.1 Cleaning

Periodically clean the weighing pan, draft shield element, bottom plate, draft shield (depending on the model), housing and terminal of your balance using a damp cloth. The maintenance interval depends on your standard operating procedure (SOP).

Please observe the following notes

WARNING



Damage of balance

- The balance must be disconnected from the power supply.
- Ensure that no liquid comes into contact with the balance, the terminal or the AC adapter.

- Never open the balance, terminal or AC adapter – they contain no components, which can be cleaned, repaired or replaced by the user.



CAUTION

Damage of balance

On no account use cleaning agents which contain solvents or abrasive ingredients, as this can result in damage to the terminal overlay.

Cleaning

Your balance is made from high quality, resistant materials and can therefore be cleaned with a commercially available, mild cleaning agent.

- 1 Remove the draft shield for thoroughly cleaning the draft shield glasses.
- 2 When you replace these parts, make sure they are in the correct position.

Note


Please contact your METTLER TOLEDO dealer for details of the available service options. Regular servicing by an authorized service engineer ensures constant accuracy for years to come and prolongs the service life of your balance.

6.2 Cleaning the Draft Shield (0.1 mg and 1 mg Models)

- 1 Remove the weighing pan, draft-shield element (0.1 mg models), pan support (1 mg models).
- 2 Lift the draft shield off the balance and place it on a clean surface.
- 3 Remove the bottom plate.
- 4 Push all the glasses back as far as they will go.
- 5 Turn the cover (1) to the front.
- 6 Pull the top glass toward the back and off.
- 7 Pull the side glasses toward the back and off. Attention

Always hold the 2 **parallel** guided glasses (side glasses and top glasses) together with one hand.


8 Clean all parts then reassemble the draft shield in the reverse order.

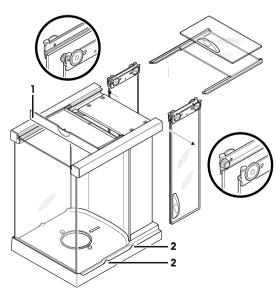

Insert glasses

Attention

Always hold the 2 **parallel** guided glasses (side glasses and top glasses) together and parallel with one hand. The side glasses must not be placed oust-side the projections (2).

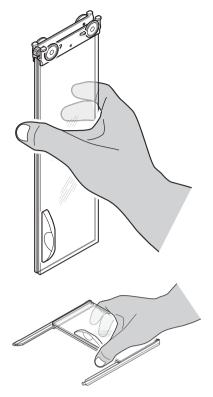

6.3 Disposal

In conformance with the European Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

Thank you for your contribution to environmental protection.



7 Technical Data

7.1 General Data "S" and "M" Weighing Platform



CAUTION

Power supply

Power supply connector with AC/DC adapter:

11107909 Primary: 10

Cable to AC adapter:

Primary: 100-240 VAC, -15%/+10%, 50/60 Hz Secondary: 12 VDC ±/-3%, 2.0 A (with electronic overload protection) Design: 3-core, with country-specific plug **Note** Make sure the power supply plug is freely accessible 12 VDC ±/-3%, 2.0 A, maximum ripple: 80 mVDCpp

Power supply to the balance:

Protection and standards

Overvoltage category: Degree of pollution: Protection:

Standards for safety and EMC: Range of application:

Environmental conditions

Height above mean sea level: Ambient temperature: Relative air humidity:

Materials

Housing: Terminal: Weighing pan: Draft shield: Draft shield element: Class II

2

IP54, in use with weighing pan inserted, protected against dust and water See Declaration of Conformity For use only in closed interior rooms

Up to 4000 m 5-40 °C Max. 80% at 31 °C, linearly decreasing to 50% at 40 °C, noncondensing

Die-cast aluminum, laquered, plastic and chrome steel Die-cast zinc, chromed and plastics Chrome-nickel steel X2CrNiMo-17-12-2 Aluminium, plastic, chrome steel and glas Die-cast zinc, chromed (10 mg models, S-platform) Chrome steel X2CrNiMo-17-13-2 (0.1 mg models)


7.2 Explanatory Notes for the METTLER TOLEDO AC Adapter

The certified external power supply which conforms to the requirements for Class II double insulated equipment is not provided with a protective earth connection but with a functional earth connection for EMC purposes. This earth connection IS NOT a safety feature. Further information about conformance of our products can be found in the brochure "Declaration of Conformity" which is coming with each product.

In case of testing with regard to the European Directive 2001/95/EC the power supply and the balance have to be handled as Class II double insulated equipment.

Consequently an earth bonding test is not required. Similarly it is not necessary to carry out an earth bonding test between the supply earth conductor and any exposed metalwork on the balance.

Because the balance are sensitive to static charges a leakage resistor, typically 10 k Ω , is connected between the earth connector and the power supply output terminals. The arrangement is shown in the equivalent circuit diagram. This resistor is not part of the electrical safety arrangement and does not require testing at regular intervals.

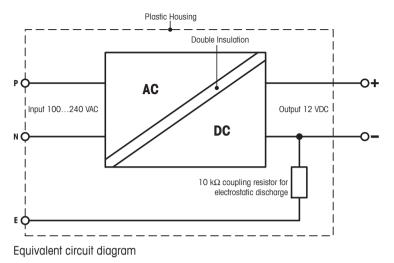

7.3 General Data "L" Weighing Platform

Power supply

Power supply:	115-240 VAC, -15%/+10%, 50/60 Hz, 0.4 A
Power cable:	3-core, with country-specific plug
Protection and standards	
Overvoltage category:	Class II
Degree of pollution:	2
Protection:	IP54, in use with weighing pan inserted, protected against dust and water
Standards for safety and EMC:	See Declaration of Conformity
Range of application:	For use only in closed interior rooms
Environmental conditions	
Environmental conditions Height above mean sea level:	Up to 4000 m
	Up to 4000 m 5-40 °C
Height above mean sea level:	
Height above mean sea level: Ambient temperature:	5-40 °C Max. 80% at 31 °C, linearly decreasing to 50% at 40 °C, non-
Height above mean sea level: Ambient temperature: Relative air humidity:	5-40 °C Max. 80% at 31 °C, linearly decreasing to 50% at 40 °C, non-
Height above mean sea level: Ambient temperature: Relative air humidity: Materials	5-40 °C Max. 80% at 31 °C, linearly decreasing to 50% at 40 °C, non- condensing
Height above mean sea level: Ambient temperature: Relative air humidity: Materials Housing:	5-40 °C Max. 80% at 31 °C, linearly decreasing to 50% at 40 °C, non- condensing Sheet aluminum, die cast, laquered, plastic and chrome steel

7.4 Model-specif Data

7.4.1 Balances with Readability of 0.1 mg, "S" Platform with Draft Shield

		XP204S	XP404S	XP404SDR
Limit values				
Maximum capacity		210 g	410 g	410 g
Readability		0.1 mg	0.1 mg	1 mg
Tare range (fromto)		0 210 g	0 410 g	0 410 g
Maximum capacity, fine range		-	_	80 g
Readability, fine range		-	_	0.1 mg
Repeatability (at nominal load)	sd	0.2 mg	0.1 mg	0.6 mg
Repeatability, fine range (at nominal load)	sd	-	_	0.1 mg
Linearity deviation		0.2 mg	0.2 mg	0.6 mg
Eccentricity deviation (test load)		0.3 mg (100 g)	0.3 mg (200 g)	1 mg (200 g)
Sensitivity offset (test weight)		1 mg (200 g)	2 mg (400 g)	2 mg (400 g)
Sensitivity temperature drift 1)		0.00015%/°C	0.00015%/°C	0.00015%/°C
Sensitivity stability		0.00025%/a	0.00025%/a	0.00025%/a
Typical values		I	1	I
Repeatability	sd	0.12 mg	0.06 mg	4 mg
Repeatability, fine range	sd	-	-	0.06 mg
Linearity deviation		0.07 mg	0.07 mg	0.07 mg
Eccentric deviation (test load)		0.08 mg (100 g)	0.08 mg (200 g)	0.08 mg (200 g)
Sensitivity offset (test load)		0.4 mg (200 g)	0.48 mg (400 g)	0.48 mg (400 g)
Minimum sample weight (according to USP)		360 mg	180 mg	1200 mg
Minimum sample weight (according to USP),		-	_	18 mg
fine range				5
Minimum sample weight (U=1%, k=2)		24 mg	12 mg	80 mg
Minimum sample weight (U=1%, k=2), fine range		_	-	12 mg
Settling time		2 s	2 s	2 s
Dimensions	1	I	1	
Balance dimensions (WxDxH)		214x395x363 mm	214x395x363 mm	214x395x363 mm
Weighing pan dimensions		90 mm (Ø)	90 mm (Ø)	90 mm (Ø)
Typical uncertainties and supplementary data				
Repeatability	sd	0.12mg+0.000015%·R- gr	0.06mg+0.000005%·R- gr	4mg+0.000025%.Rgr
Repeatability, fine range	sd	_	_	0.06mg+0.000025%·R gr
Differential linearity deviation	sd	√(6pg·Rnt)	√(3pg·Rnt)	√(3pg·Rnt)
Differential eccentric load deviation	sd	0.00004%·Rnt	0.00002%·Rnt	0.00002%·Rnt
Sensitivity offset	sd	0.0001%·Rnt	0.00006%·Rnt	0.00006%·Rnt
Minimum sample weight (according to USP)		360mg+0.045%.Rgr	180mg+0.015%.Rgr	1200mg+0.075%.Rgr
Minimum sample weight (according to USP), fine range		_	-	18mg+0.075%-Rgr
Minimum sample weight (U=1%, k=2)		24mg+0.003%.Rgr	12mg+0.001%.Rgr	80mg+0.005%·Rgr
Minimum sample weight (U=1%, k=2), fine range		_	-	12mg+0.005%·Rgr
Interface update rate		23 /s	23 /s	23 /s
Usable height of draft shield		248 mm	248 mm	248 mm
Weight of balance		8.2 kg	8.2 kg	8.2 kg
Number of built-in reference weights	1	1	1	1
Weights for routine testing	1	1	1	1
OIML CarePac		200 g F2, 10 g F1	200 g F2, 20 g F1	200 g F2, 20 g F1
	Voiabta			
	veignis	#11123001	#11123000	#11123000
ASTM CarePac		200 g 1, 10 g 1	200 g 1, 20 g 1	200 g 1, 20 g 1
,		#11123101	#11123100	#11123100

sd = Standard deviation

Rgr = Gross weight

- ¹⁾ In the temperature range 10...30 °C
- Rnt = Net weight (sample weight)
 - = Year (annum)

α

7.4.2 Balances with Readability of 1 mg, "S" Platform with Draft Shield

		XP203S	XP603S	XP603SDR
Limit values			-	
Maximum capacity		210 g	610 mg	610 g
Readability		1 mg	1 mg	10 mg
Tare range (fromto)		0 210 g	0 610 g	0 610 g
Maximum capacity, fine range		_	-	120 g
Readability, fine range		-	-	1 mg
Repeatability (at nominal load)	sd	0.9 mg	0.9 mg	6 mg
Repeatability, fine range (at nominal load)	sd	-	-	1 mg
Linearity deviation		2 mg	2 mg	6 mg
Eccentricity deviation (test load)		3 mg (200 g)	3 mg (200 g)	10 mg (200 g)
Sensitivity offset (test weight)		5 mg (200 g)	4.5 mg (600 g)	9 mg (600 g)
Sensitivity temperature drift 1)		0.0005%/°C	0.0002%/°C	0.0002%/°C
Sensitivity stability		0.0025%/a	0.001%/a	0.001%/a
Typical values			-	
Repeatability	sd	0.5 mg	0.5 mg	4 mg
Repeatability, fine range	sd	_	-	0.8 mg
Linearity deviation		0.65 mg	0.7 mg	0.7 mg
Eccentric deviation (test load)		0.6 mg (200 g)	0.8 mg (200 g)	0.8 mg (200 g)
Sensitivity offset (test load)		3.2 mg (400 g)	2.4 mg (600 g)	6 mg (600 g)
Minimum sample weight (according to USP)		1500 mg	1.5 g	12 g
Minimum sample weight (according to USP), ine range		-	-	2.4 g
Vinimum sample weight (U=1%, k=2)		100 mg	100 mg	800 mg
Vinimum sample weight (U=1%, k=2), fine range		_	-	160 mg
Settling time		1.5 s	1.5 s	1.5 s
Dimensions	1			
Balance dimensions (WxDxH)		198x392x276 mm	198x392x276 mm	198x392x276 mm
Weighing pan dimensions		127x127 mm (WxD)	127x127 mm (WxD)	127x127 mm (WxD)
Typical uncertainties and supplementary data				
Repeatability	sd	0.5mg+0.00008%.Rgr	0.5mg+0.000025%·Rgr	4mg+0.00015%.Rgr
Repeatability, fine range	sd	_	_	0.8mg
Differential linearity deviation	sd	√(500pg⋅Rnt)	√(200pg·Rnt)	$\sqrt{(200 \text{pg} \cdot \text{Rnt})}$
Differential eccentric load deviation	sd	0.00015%·Rnt	0.0002%·Rnt	0.0002%·Rnt
Sensitivity offset	sd	0.0008%·Rnt	0.0002%·Rnt	0.0005%·Rnt
Minimum sample weight (according to USP)	1	1500mg+0.24%·Rgr	1.5g+0.075%·Rgr	12g+0.45%·Rgr
Minimum sample weight (according to USP), ine range		_	_	2.4g
Minimum sample weight (U=1%, k=2)		100mg+0.016%.Rgr	100mg+0.005%·Rgr	800mg+0.03%·Rgr
Minimum sample weight (U=1%, k=2), fine range		_	-	160mg
Interface update rate	1	23 /s	23 /s	23 /s
Jsable height of draft shield		175 mm	175 mm	175 mm
Weight of balance		7.7 kg	7.7 kg	7.7 kg
Number of built-in reference weights		1	1	1
	1	1		1
Weights for routine testing DIML CarePac		200 g F2, 10 g F1	500 g F2, 20 g F1	500 g F2, 20 g F1

	XP203S	XP603S	XP603SDR
ASTM CarePac	200 g 1, 10 g 1	500 g 1, 20 g 1	500 g 1, 20 g 1
Weights	#11123101	#11123107	#11123107

α

sd = Standard deviation

Rgr = Gross weight

Rnt = Net weight (sample weight)

= Year (annum)

¹⁾ In the temperature range 10...30 °C

		XP1203S	XP2003SDR	XP5003SDR
Limit values				
Maximum capacity		1210 g	2.1 kg	5.1 kg
Readability		1 mg	10 mg	10 mg
Tare range (fromto)		0 1210 g	0 2.1 kg	0 5.1 kg
Maximum capacity, fine range		-	0.5 kg	1 kg
Readability, fine range		-	1 mg	1 mg
Repeatability (at nominal load)	sd	0.8 mg	6 mg	6 mg
Repeatability, fine range (at nominal load)	sd	-	1 mg	1 mg
Linearity deviation		2 mg	6 mg	6 mg
Eccentricity deviation (test load)		3 mg (500 g)	10 mg (1 kg)	10 mg (2 kg)
Sensitivity offset (test weight)		6 mg (1200 g)	10 mg (2 kg)	20 mg (5 kg)
Sensitivity temperature drift 1)		0.0002%/°C	0.0003%/°C	0.0003%/°C
Sensitivity stability		0.001%/a	0.0025%/a	0.0015%/a
Typical values	-	1	1	1
Repeatability	sd	0.4 mg	4 mg	4 mg
Repeatability, fine range	sd	-	0.6 mg	0.6 mg
Linearity deviation	1	0.7 mg	0.7 mg	1 mg
Eccentric deviation (test load)		1 mg (500 g)	0.6 mg (1 kg)	0.6 mg (2 kg)
Sensitivity offset (test load)		2.9 mg (1200 g)	3.2 mg (2 kg)	10 mg (5 kg)
Minimum sample weight (according to USP)		1.2 g	12 g	12 g
Minimum sample weight (according to USP), fine range		-	1.8 g	1.8 g
Minimum sample weight (U=1%, k=2)		80 mg	800 mg	800 mg
Minimum sample weight (U=1%, k=2), fine range		-	120 mg	120 mg
Settling time		1.5 s	2 s	2 s
Dimensions				
Balance dimensions (WxDxH)		198x392x276 mm	214x395x363 mm	214x395x363 mm
Weighing pan dimensions		127x127 mm (WxD)	127x127 mm (WxD)	127x127 mm (WxD)
Typical uncertainties and supplementary data				
Repeatability	sd	0.4mg+0.000015%.Rgr	4mg+0.00005%.Rgr	4mg+0.00002%.Rgr
Repeatability, fine range	sd	-	0.6mg+0.00004%·Rgr	0.6mg+0.00002%·Rgr
Differential linearity deviation	sd	√(100pg·Rnt)	√(60pg·Rnt)	√(50pg·Rnt)
Differential eccentric load deviation	sd	0.0001%·Rnt	0.00003%·Rnt	0.000015%·Rnt
Sensitivity offset	sd	0.00012%·Rnt	0.00008%·Rnt	0.0001%·Rnt
Minimum sample weight (according to USP)		1.2g+0.05%·Rgr	12g+0.15%·Rgr	12g+0.06%·Rgr
Minimum sample weight (according to USP), fine range		-	1.8g+0.12%·Rgr	1.8g+0.06%.Rgr
Minimum sample weight (U=1%, k=2)		80mg+0.003%.Rgr	800mg+0.01%.Rgr	800mg+0.004%·Rgr
Minimum sample weight (U=1%, k=2), fine range		-	120mg+0.008%.Rgr	120mg+0.004%·Rgr
Interface update rate		23 /s	23 /s	23 /s
Usable height of draft shield		175 mm	-	-
Weight of balance		7.7 kg	8.6 kg	8.6 kg
Number of built-in reference weights		1	1	1
Weights for routine testing				
OIML CarePac		1000 g F2, 50 g F2	2000 g F2, 100 g F2	5000 g F2, 200 g F2
		-	-	-

				XP1203S		XP2003SDR	XP5003SDR
ASTN	1 CareF	Pac		1000 g 1, 50 g 1		2000 g 1, 100 g 1	5000 g 1, 200 g 1
			Weights	#11123108		#11123109	#11123111
sd	=	Standard deviation		Rnt	=	Net weight (sample weight)	
Rgr	=	Gross weight		a	=	Year (annum)	

¹⁾ In the temperature range 10...30 °C

7.4.3 Balances with Readability of 10 mg, "S" Platform with Draft Shield Element

		XP802S	XP1202S	XP2002S
Limit values				
Maximum capacity		810 g	1210 g	2.1 kg
Readability		10 mg	10 mg	10 mg
Tare range (fromto)		0 810 g	0 1210 g	0 2.1 kg
Repeatability (at nominal load)	sd	8 mg	8 mg	8 mg
Linearity deviation		20 mg	20 mg	20 mg
Eccentricity deviation (test load)		20 mg (500 g)	20 mg (500 g)	30 mg (1 kg)
Sensitivity offset (test weight)		60 mg (800 g)	60 mg (1200 g)	60 mg (2 kg)
Sensitivity temperature drift 1)		0.0003%/°C	0.0003%/°C	0.0003%/°C
Sensitivity stability		0.0025%/a	0.0025%/a	0.0025%/a
Typical values		•		·
Repeatability	sd	4 mg	4 mg	4 mg
Linearity deviation		7 mg	7 mg	7 mg
Eccentric deviation (test load)		3 mg (500 g)	3 mg (500 g)	3 mg (1 kg)
Sensitivity offset (test load)		32 mg (800 g)	24 mg (1200 g)	32 mg (2 kg)
Minimum sample weight (according to USP)		12 g	12 g	12 g
Minimum sample weight (U=1%, k=2)		800 mg	800 mg	800 mg
Settling time		1.2 s	1.2 s	1.2 s
Dimensions			1	
Balance dimensions (WxDxH)		194x392x96 mm	194x392x96 mm	194x392x96 mm
Weighing pan dimensions		170x205 mm (WxD)	170x205 mm (WxD)	170x205 mm (WxD)
Typical uncertainties and supplementary data		·		·
Repeatability	sd	4mg+0.00025%.Rgr	4mg+0.00015%.Rgr	4mg+0.0001%.Rgr
Differential linearity deviation	sd	√(15ng·Rnt)	√(10ng⋅Rnt)	√(6ng⋅Rnt)
Differential eccentric load deviation	sd	0.0003%·Rnt	0.0003%·Rnt	0.00015%·Rnt
Sensitivity offset	sd	0.002%·Rnt	0.001%·Rnt	0.0008%·Rnt
Minimum sample weight (according to USP)		12g+0.75%·Rgr	12g+0.45%·Rgr	12g+0.3%·Rgr
Minimum sample weight (U=1%, k=2)		800mg+0.05%·Rgr	800mg+0.03%.Rgr	800mg+0.02%.Rgr
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		7.1 kg	6.6 kg	7.1 kg
Number of built-in reference weights		1	1	1
Weights for routine testing	-			
OIML CarePac		500 g F2, 20 g F1	1000 g F2, 50 g F2	2000 g F2, 100 g F2
V	Veights	#11123007	#11123008	#11123009
ASTM CarePac		500 g 1, 20 g 1	1000 g 1, 50 g 1	2000 g 1, 100 g 1
	Voiabto	#11123107	#11123108	#11123109

Rgr = Gross weight

a = Year (annum)

¹⁾ In the temperature range 10...30 °C

	XP4002S	XP4002SDR	XP6002S
Limit values			
Maximum capacity	4.1 kg	4.1 kg	6.1 kg
Readability	10 mg	100 mg	10 mg

		XP4002S	XP4002SDR	XP6002S
Tare range (fromto)		0 4.1 kg	0 4.1 kg	0 6.1 kg
Maximum capacity, fine range		-	0.8 kg	-
Readability, fine range		-	10 mg	-
Repeatability (at nominal load)	sd	8 mg	80 mg	8 mg
Repeatability, fine range (at nominal load)	sd	_	8 mg	-
Linearity deviation		20 mg	60 mg	20 mg
Eccentricity deviation (test load)		30 mg (2 kg)	100 mg (2 kg)	30 mg (2 kg)
Sensitivity offset (test weight)		60 mg (4 kg)	60 mg (4 kg)	60 mg (6 kg)
Sensitivity temperature drift 1)		0.0003%/°C	0.0003%/°C	0.0003%/°C
Sensitivity stability		0.0015%/a	0.0015%/a	0.0015%/a
Typical values				
Repeatability	sd	4 mg	40 mg	4 mg
Repeatability, fine range	sd	-	40 mg	-
Linearity deviation		7 mg	7 mg	7 mg
Eccentric deviation (test load)		8 mg (2 kg)	6 mg (2 kg)	8 mg (2 kg)
Sensitivity offset (test load)		32 mg (4 kg)	32 mg (4 kg)	30 mg (6 kg)
Minimum sample weight (according to USP)		12 g	120 g	12 g
Minimum sample weight (according to USP), fine range		_	12 g	_
Minimum sample weight (U=1%, k=2)		800 mg	8 g	800 mg
Minimum sample weight (U=1%, k=2), fine range		-	0.8 g	-
Settling time		1.2 s	1.2 s	1.2 s
Dimensions				•
Balance dimensions (WxDxH)		194x392x96 mm	194x392x96 mm	194x392x96 mm
Weighing pan dimensions		170x205 mm (WxD)	170x205 mm (WxD)	170x205 mm (WxD)
Typical uncertainties and supplementary dat	a	-		
Repeatability	sd	4mg+0.00005%·Rgr	40mg+0.00025%.Rgr	4mg+0.00003%.Rgr
Repeatability, fine range	sd	-	40mg+0.00025%.Rgr	-
Differential linearity deviation	sd	√(3ng⋅Rnt)	√(3ng⋅Rnt)	√(2ng⋅Rnt)
Differential eccentric load deviation	sd	0.0002%·Rnt	0.00015%·Rnt	0.0002%·Rnt
Sensitivity offset	sd	0.0004%·Rnt	0.0004%·Rnt	0.00025%·Rnt
Minimum sample weight (according to USP)		12g+0.15%·Rgr	120g+0.75%·Rgr	12g+0.09%·Rgr
Minimum sample weight (according to USP), fine range		_	12g+0.75%.Rgr	-
Minimum sample weight (U=1%, k=2)		800mg+0.01%·Rgr	8g+0.05%·Rgr	800mg+0.006%·Rgr
Minimum sample weight (U=1%, k=2), fine range		_	0.8g+0.05%·Rgr	_
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		6.6 kg	7.1 kg	6.6 kg
Number of built-in reference weights		1	1	1
Weights for routine testing				
OIML CarePac		2000 g F2, 200 g F2	2000 g F2, 200 g F2	5000 g F2, 200 g F2
	Weights	#11123010	#11123010	#11123011
ASTM CarePac		2000 g 4, 200 g 4	2000 g 4, 200 g 4	5000 g 4, 200 g 4
	Weights	#11123110	#11123110	#11123111
sd = Standard deviation		Rnt =	Net weight (sample weigh	t)
Rgr = Gross weight		a =	Year (annum)	

 $^{1)}$ In the temperature range 10...30 $^{\circ}\text{C}$

	XP6002SDR	XP8002S
Limit values		
Maximum capacity	6.1 kg	8.1 kg
Readability	100 mg	10 mg
Tare range (fromto)	0 6.1 kg	0 8.1 kg

		XP6002SDR	XP8002S
Maximum capacity, fine range		1.2 kg	-
Readability, fine range		10 mg	-
Repeatability (at nominal load)	sd	60 mg	8 mg
Repeatability, fine range (at nominal load)	sd	8 mg	-
Linearity deviation		60 mg	20 mg
Eccentricity deviation (test load)		100 mg (2 kg)	40 mg (5 kg)
Sensitivity offset (test weight)		150 mg (6 kg)	60 mg (8 kg)
Sensitivity temperature drift 1)		0.0003%/°C	0.00025%/°C
Sensitivity stability		0.0015%/a	0.0015%/a
Typical values		1	-
Repeatability	sd	40 mg	4 mg
Repeatability, fine range	sd	5 mg	-
Linearity deviation		7 mg	7 mg
Eccentric deviation (test load)		8 mg (2 kg)	10 mg (5 kg)
Sensitivity offset (test load)		30 mg (6 kg)	32 mg (8 kg)
Minimum sample weight (according to USP)	1	120 g	12 g
Minimum sample weight (according to USP),		15 g	-
fine range			
Minimum sample weight (U=1%, k=2)		8 g	800 mg
Minimum sample weight		1 g	-
(U=1%, k=2), fine range			
Settling time		1.2 s	1.5 s
Dimensions	-	1	1
Balance dimensions (WxDxH)		194x392x96 mm	194x392x96 mm
Weighing pan dimensions		170x205 mm (WxD)	170x205 mm (WxD)
Typical uncertainties and supplementary data			
Repeatability	sd	40mg+0.00015%·Rgr	4mg+0.000025%·Rg
Repeatability, fine range	sd	5mg	-
Differential linearity deviation	sd	√(2ng⋅Rnt)	√(1.5ng·Rnt)
Differential eccentric load deviation	sd	0.0002%·Rnt	0.0001%·Rnt
Sensitivity offset	sd	0.00025%·Rnt	0.0002%·Rnt
Minimum sample weight (according to USP)		120g+0.45%·Rgr	12g+0.075%.Rgr
Minimum sample weight (according to USP), fine range		15g	-
Minimum sample weight (U=1%, k=2)		8g+0.03%·Rgr	800mg+0.005%·Rgr
Minimum sample weight (U=1%, k=2), fine range		1g	-
Interface update rate		23 /s	23 /s
Weight of balance		6.6 kg	6.6 kg
Number of built-in reference weights		1	1
Weights for routine testing			
OIML CarePac		5000 g F2, 200 g F2	5000 g F2, 200 g F2
V	Veights	#11123011	#11123011
ASTM CarePac		5000 g 4, 200 g 4	5000 g 4, 200 g 4
	Veiahts	#11123111	#11123111
sd = Standard deviation Rgr = Gross weight		Rnt = a =	Net weight (sample wei Year (annum)

¹⁾ In the temperature range 10...30 °C

	XP10002S	XP10002SDR	XP12002SDR		
Limit values					
Maximum capacity	10.1 kg	10.1 kg	12.1 kg		
Readability	10 mg	100 mg	100 mg		
Tare range (fromto)	0 10.1 kg	0 10.1 kg	0 12.1 kg		
Maximum capacity, fine range	-	2 kg	2.4 kg		

		XP10002S	XP10002SDR	XP12002SDR
Readability, fine range		-	10 mg	10 mg
Repeatability (at nominal load)	sd	8 mg	60 mg	60 mg
Repeatability, fine range (at nominal load)	sd	-	8 mg	10 mg
Linearity deviation		20 mg	50 mg	60 mg
Eccentricity deviation (test load)		40 mg (5 kg)	100 mg (5 kg)	100 mg (5 kg)
Sensitivity offset (test weight)		50 mg (10 kg)	100 mg (10 kg)	96 mg (12 kg)
Sensitivity temperature drift 1)		0.00025%/°C	0.00025%/°C	0.00025%/°C
Sensitivity stability		0.0015%/a	0.0015%/a	0.00015%/a
Typical values				·
Repeatability	sd	4 mg	40 mg	40 mg
Repeatability, fine range	sd	-	5 mg	6 mg
Linearity deviation		6.5 mg	4 mg	7 mg
Eccentric deviation (test load)		10 mg (5 kg)	10 mg (5 kg)	10 mg (5 kg)
Sensitivity offset (test load)		30 mg (10 kg)	30 mg (10 kg)	60 mg (12 kg)
Minimum sample weight (according to USP)	1	12 g	120 g	120 g
Minimum sample weight (according to USP), fine range		_	15 g	18 g
Minimum sample weight (U=1%, k=2)		800 mg	8 g	8 g
Minimum sample weight (U=1%, k=2), fine range		-	1 g	1.2 g
Settling time		1.5 s	1.5 s	1.8 s
Dimensions			L	
Balance dimensions (WxDxH)		194x392x96 mm	194x392x96 mm	194x392x96 mm
Weighing pan dimensions		170x205 mm (WxD)	170x205 mm (WxD)	170x205 mm (WxD)
Typical uncertainties and supplementary date	1	1		
Repeatability	sd	4mg+0.00002%.Rgr	40mg+0.0001%.Rgr	40mg+0.00008%.Rg
Repeatability, fine range	sd	-	5mg	6mg+0.00008%.Rgr
Differential linearity deviation	sd	√(1ng·Rnt)	√(400pg⋅Rnt)	√(1ng·Rnt)
Differential eccentric load deviation	sd	0.0001%·Rnt	0.0001%·Rnt	0.0001%·Rnt
Sensitivity offset	sd	0.00015%·Rnt	0.00015%·Rnt	0.00025%·Rnt
Minimum sample weight (according to USP)		12g+0.06%·Rgr	120g+0.3%·Rgr	120g+0.24%·Rgr
Minimum sample weight (according to USP), fine range		-	15g	18g+0.24%·Rgr
Minimum sample weight (U=1%, k=2)		800mg+0.004%·Rgr	8g+0.02%.Rgr	8g+0.016%·Rgr
Minimum sample weight (U=1%, k=2), fine range		-	1g	1.2g+0.016%·Rgr
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		6.6 kg	6.6 kg	8.1 kg
Number of built-in reference weights		1	1	1
Weights for routine testing				
OIML Single Weights		500 g F2	500 g F2	500 g F2
	Small	#11118203	#11118203	#11118203
OIML Single Weights		10000 g F2	10000 g F2	10000 g F2
	Large	#11118211	#11118211	#11118211
sd = Standard deviation		Rnt =	Net weight (sample weig	nt)
Rgr = Gross weight		a =	Year (annum)	

¹⁾ In the temperature range 10...30 °C

7.4.4 Balances with Readability of 0.1 g, "S" Platform

	XP2001S	XP4001S	XP6001S		
Limit values					
Maximum capacity	2.1 kg	4.1 kg	6.1 kg		
Readability	100 mg	100 mg	100 mg		
Tare range (fromto)	0 2.1 kg	0 4.1 kg	0 6.1 kg		

		XP2001S	XP4001S	XP6001S
Repeatability (at nominal load)	sd	80 mg	80 mg	80 mg
Linearity deviation		60 mg	60 mg	60 mg
Eccentricity deviation (test load)		100 mg (1 kg)	200 mg (2 kg)	200 mg (2 kg)
Sensitivity offset (test weight)		150 mg (2 kg)	240 mg (4 kg)	240 mg (6 kg)
Sensitivity temperature drift 1)		0.0015%/°C	0.0015%/°C	0.0015%/°C
Sensitivity stability		0.005%/a	0.005%/a	0.005%/a
Typical values			L	·
Repeatability	sd	40 mg	40 mg	40 mg
Linearity deviation		20 mg	13 mg	19 mg
Eccentric deviation (test load)		20 mg (1 kg)	32 mg (2 kg)	32 mg (2 kg)
Sensitivity offset (test load)		80 mg (2 kg)	160 mg (4 kg)	140 mg (6 kg)
Minimum sample weight (according to USP)		120 g	120 g	120 g
Minimum sample weight (U=1%, k=2)		8 g	8 g	8 g
Settling time		0.8 s	0.8 s	0.8 s
Dimensions			·	
Balance dimensions (WxDxH)		194x392x96 mm	194x392x96 mm	194x392x96 mm
Weighing pan dimensions		190x223 mm (WxD)	190x223 mm (WxD)	190x223 mm (WxD)
Typical uncertainties and supplementary data	1			
Repeatability	sd	40mg+0.001%·Rgr	40mg+0.0005%.Rgr	40mg+0.0003%.Rgr
Differential linearity deviation	sd	√(50ng·Rnt)	√(10ng·Rnt)	√(15ng·Rnt)
Differential eccentric load deviation	sd	0.001%·Rnt	0.0008%·Rnt	0.0008%·Rnt
Sensitivity offset	sd	0.002%·Rnt	0.002%·Rnt	0.0012%·Rnt
Minimum sample weight (according to USP)		120g+3%·Rgr	120g+1.5%·Rgr	120g+0.9%.Rgr
Minimum sample weight (U=1%, k=2)		8g+0.2%·Rgr	8g+0.1%·Rgr	8g+0.06%·Rgr
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		6.6 kg	6.6 kg	6.6 kg
Number of built-in reference weights		1	1	1
Weights for routine testing				
OIML CarePac		2000 g F2, 100 g F2	2000 g F2, 200 g F2	5000 g F2, 200 g F2
	Weights	#11123009	#11123010	#11123011
ASTM CarePac		2000 g 1, 100 g 1	2000 g 4, 200 g 4	5000 g 4, 200 g 4
		#11123109	#11123110	#11123111

α

Standard deviation sd =

Year (annum) =

Gross weight Rgr =

1) In the temperature range 10...30 $^\circ\text{C}$

		XP8001S	XP10001S
Limit values			
Maximum capacity		8.1 kg	10.1 kg
Readability		100 mg	100 mg
Tare range (fromto)		0 8.1 kg	0 10.1 kg
Repeatability (at nominal load)	sd	80 mg	80 mg
Linearity deviation		100 mg	100 mg
Eccentricity deviation (test load)		200 mg (5 kg)	200 mg (5 kg)
Sensitivity offset (test weight)		600 mg (8 kg)	500 mg (10 kg)
Sensitivity temperature drift 1)		0.0015%/°C	0.0015%/°C
Sensitivity stability		0.005%/a	0.005%/a
Typical values			
Repeatability	sd	40 mg	40 mg
Linearity deviation		34 mg	34 mg
Eccentric deviation (test load)		30 mg (5 kg)	30 mg (5 kg)
Sensitivity offset (test load)		320 mg (8 kg)	300 mg (10 kg)
Minimum sample weight (according to USP)		120 g	120 g
Minimum sample weight (U=1%, k=2)		8 g	8 g

		XP8001S	XP10001S
Settling time		1 s	1 s
Dimensions			
Balance dimensions (WxDxH)		194x392x96 mm	194x392x96 mm
Weighing pan dimensions		190x223 mm (WxD)	190x223 mm (WxD)
Typical uncertainties and supplementary date	1		
Repeatability	sd	40mg+0.00025%·Rgr	40mg+0.0002%.Rgr
Differential linearity deviation	sd	√(35ng·Rnt)	√(30ng·Rnt)
Differential eccentric load deviation	sd	0.0003%·Rnt	0.0003%·Rnt
Sensitivity offset	sd	0.002%·Rnt	0.0015%·Rnt
Minimum sample weight (according to USP)		120g+0.75%.Rgr	120g+0.6%·Rgr
Minimum sample weight (U=1%, k=2)		8g+0.05%∙Rgr	8g+0.04%·Rgr
Interface update rate		23 /s	23 /s
Weight of balance		6.6 kg	6.6 kg
Number of built-in reference weights		1	1
Weights for routine testing			
OIML CarePac		5000 g F2, 200 g F2	500 g F2
, in the second s	Weights	#11123011	#11118203
			OIML Single Weights, Small
ASTM CarePac		5000 g 4, 200 g 4	10000 g F2
l l l l l l l l l l l l l l l l l l l	Weights	#11123111	#11118211
			OIML Single Weights, Large
sd = Standard deviation		Rnt =	Net weight (sample weight
Rgr = Gross weight		a =	Year (annum)

¹⁾ In the temperature range 10...30 °C

7.4.5 Balances with Readability of 10 mg / 0.1 g / 1 g, "M" Platform

		XP6002MDR	XP12002MDR	XP8001M
Limit values		1		I
Maximum capacity		6.1 kg	12.1 kg	8.1 kg
Readability		100 mg	100 mg	100 mg
Tare range (fromto)		0 6.1 kg	0 12.1 kg	0 8.1 kg
Maximum capacity, fine range		1.2 kg	2.4 kg	-
Readability, fine range		10 mg	10 mg	-
Repeatability (at nominal load)	sd	60 mg	60 mg	80 mg
Repeatability, fine range (at nominal load)	sd	10 mg	10 mg	-
Linearity deviation		60 mg	60 mg	100 mg
Eccentricity deviation (test load)		100 mg (2 kg)	100 mg (5 kg)	200 mg (5 kg)
Sensitivity offset (test weight)		150 mg (6 kg)	96 mg (12 kg)	600 mg (8 kg)
Sensitivity temperature drift 1)		0.0003%/°C	0.00025%/°C	0.0015%/°C
Sensitivity stability		0.0015%/a	0.0015%/a	0.005%/a
Typical values				
Repeatability	sd	40 mg	40 mg	40 mg
Repeatability, fine range	sd	6 mg	6 mg	-
Linearity deviation		7 mg	7 mg	36 mg
Eccentric deviation (test load)		10 mg (2 kg)	10 mg (5 kg)	30 mg (5 kg)
Sensitivity offset (test load)		60 mg (6 kg)	60 mg (12 kg)	320 mg (8 kg)
Minimum sample weight (according to USP)		120 g	120 g	120 g
Minimum sample weight (according to USP), fine range		18 g	18 g	-
Minimum sample weight (U=1%, k=2)		8 g	8 g	8 g

		XP6002MDR	XP12002MDR	XP8001M
Minimum sample weight (U=1%, k=2), fine range		1.2 g	1.2 g	-
Settling time		1.5 s	1.8 s	1.2 s
Dimensions				- .
Balance dimensions (WxDxH)		240x419x110 mm	240x419x110 mm	240x419x110 mm
Weighing pan dimensions		237x237 mm (WxD)	237x237 mm (WxD)	237x237 mm (WxD)
Typical uncertainties and supplementary data	I		·	
Repeatability	sd	40mg+0.00015%.Rgr	40mg+0.00008%.Rgr	40mg+0.00025%.Rgr
Repeatability, fine range	sd	6mg+0.00015%.Rgr	6mg+0.00008%.Rgr	-
Differential linearity deviation	sd	√(2ng·Rnt)	√(1ng·Rnt)	√(40ng·Rnt)
Differential eccentric load deviation	sd	0.00025%·Rnt	0.0001%·Rnt	0.0003%·Rnt
Sensitivity offset	sd	0.0005%·Rnt	0.00025%·Rnt	0.002%·Rnt
Minimum sample weight (according to USP)		120g+0.45%·Rgr	120g+0.24%.Rgr	120g+0.75%.Rgr
Minimum sample weight (according to USP), fine range		18g+0.45%·Rgr	18g+0.24%·Rgr	-
Minimum sample weight (U=1%, k=2)		8g+0.03%·Rgr	8g+0.016%.Rgr	8g+0.05%·Rgr
Minimum sample weight (U=1%, k=2), fine range		1.2g+0.03%·Rgr	1.2g+0.016%·Rgr	-
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		8.1 kg	8.1 kg	8.1 kg
Number of built-in reference weights		1	1	1
Weights for routine testing				·
OIML Single Weights		200 g F2	500 g F2	200 g F2
	Small	#11118202	#11118203	#11118202
OIML Single Weights		5000 g F2	10000 g F2	5000 g F2
	Large	#11118206	#11118211	#11118206
sd = Standard deviation		Rnt =	Net weight (sample weigh	it)
sd = Standard deviation		Rnt =	Net weight (sample weigh	it)

Rgr = Gross weight

= Year (annum)

¹⁾ In the temperature range 10...30 °C

		XP8001MDR	XP12001M	XP12000M*
Limit values				
Maximum capacity		8.1 kg	12.1 kg	12.1 kg
Readability		1000 mg	100 mg	1000 mg
Tare range (fromto)		0 8.1 kg	0 12.1 kg	0 12.1 kg
Maximum capacity, fine range		1.6 kg	-	-
Readability, fine range		100 mg	-	-
Repeatability (at nominal load)	sd	600 mg	80 mg	600 mg
Repeatability, fine range (at nominal load)	sd	80 mg	-	-
Linearity deviation		600 mg	100 mg	600 mg
Eccentricity deviation (test load)		1000 mg (5 kg)	200 mg (5 kg)	1000 mg (5 kg)
Sensitivity offset (test weight)		600 mg (8 kg)	600 mg (12 kg)	600 mg (12 kg)
Sensitivity temperature drift 1)		0.0015%/°C	0.0015%/°C	0.0015%/°C
Sensitivity stability		0.005%/a	0.005%/a	0.005%/a
Typical values			· ·	
Repeatability	sd	400 mg	40 mg	400 mg
Repeatability, fine range	sd	40 mg	-	-
Linearity deviation		36 mg	34 mg	34 mg
Eccentric deviation (test load)		30 mg (5 kg)	30 mg (5 kg)	30 mg (5 kg)
Sensitivity offset (test load)		320 mg (8 kg)	290 mg (12 kg)	290 mg (12 kg)
Minimum sample weight (according to USP)		1200 g	120 g	1200 g
Minimum sample weight (according to USP), fine range		120 g	-	-
Minimum sample weight (U=1%, k=2)		80 g	8 g	80 g

α

		XP8001MDR	XP12001M	XP12000M*
Minimum sample weight		8 g	-	-
(U=1%, k=2), fine range				
Settling time		1.2 s	1.2 s	1 s
Dimensions				
Balance dimensions (WxDxH)		240x419x110 mm	240x419x110 mm	240x419x110 mm
Weighing pan dimensions		237x237 mm (WxD)	237x237 mm (WxD)	237x237 mm (WxD)
Typical uncertainties and supplementary date	l			
Repeatability	sd	400mg+0.0012%.Rgr	40mg+0.00015%.Rgr	400mg+0.0008%.Rgr
Repeatability, fine range	sd	40mg+0.0012%·Rgr	_	-
Differential linearity deviation	sd	√(40ng⋅Rnt)	√(25ng·Rnt)	√(25ng·Rnt)
Differential eccentric load deviation	sd	0.0003%·Rnt	0.0003%·Rnt	0.0003%·Rnt
Sensitivity offset	sd	0.002%·Rnt	0.0012%·Rnt	0.0012%·Rnt
Minimum sample weight (according to USP)		1200g+3.6%·Rgr	120g+0.45%·Rgr	1200g+2.4%·Rgr
Minimum sample weight (according to USP),		120g+3.6%·Rgr	-	-
fine range				
Minimum sample weight (U=1%, k=2)		80g+0.24%·Rgr	8g+0.03%·Rgr	80g+0.16%·Rgr
Minimum sample weight (U=1%, k=2), fine range		8g+0.24%·Rgr	-	_
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		8.1 kg	8.1 kg	8.1 kg
Number of built-in reference weights		1	1	1
Weights for routine testing			•	1
OIML Single Weights		200 g F2	500 g F2	500 g F2
	Small	#11118202	#11118203	#11118203
OIML Single Weights		5000 g F2	10000 g F2	10000 g F2
	Large	#11118206	#11118211	#11118211
sd = Standard deviation		Rnt =	Net weight (sample weigh	t)
Rgr = Gross weight		a =	Year (annum)	

1) In the temperature range 10...30 °C

		XP16001M	XP16001MDR	XP20001M
Limit values			L.	
Maximum capacity		16.1 kg	16.1 kg	20.1 kg
Readability		100 mg	1000 mg	100 mg
Tare range (fromto)		0 16.1 kg	0 16.1 kg	0 20.1 kg
Maximum capacity, fine range		-	3.2 kg	-
Readability, fine range		-	100 mg	-
Repeatability (at nominal load)	sd	80 mg	600 mg	80 mg
Repeatability, fine range (at nominal load)	sd	-	80 mg	-
Linearity deviation		200 mg	600 mg	200 mg
Eccentricity deviation (test load)		200 mg (5 kg)	1000 mg (5 kg)	200 mg (10 kg)
Sensitivity offset (test weight)		800 mg (16 kg)	800 mg (16 kg)	800 mg (20 kg)
Sensitivity temperature drift 1)		0.0015%/°C	0.0015%/°C	0.0015%/°C
Sensitivity stability		0.005%/a	0.005%/a	0.005%/a
Typical values			·	· ·
Repeatability	sd	40 mg	400 mg	40 mg
Repeatability, fine range	sd	-	40 mg	_
Linearity deviation		130 mg	130 mg	130 mg
Eccentric deviation (test load)		120 mg (5 kg)	30 mg (5 kg)	120 mg (10 kg)
Sensitivity offset (test load)		260 mg (16 kg)	260 mg (16 kg)	240 mg (20 kg)
Minimum sample weight (according to USP)		120 g	1200 g	120 g
Minimum sample weight (according to USP), fine range		_	120 g	-
Minimum sample weight (U=1%, k=2)		8 g	80 g	8 g

		XP16001M	XP16001MDR	XP20001M
Minimum sample weight (U=1%, k=2), fine range		_	8 h	-
Settling time		1.2 s	1 s	1.2 s
Dimensions	1			
Balance dimensions (WxDxH)		240x419x110 mm	240x419x110 mm	240x419x110 mm
Weighing pan dimensions		237x237 mm (WxD)	237x237 mm (WxD)	237x237 mm (WxD)
Typical uncertainties and supplementary data				
Repeatability	sd	40mg+0.00012%.Rgr	400mg+0.0006%.Rgr	40mg+0.0001%.Rgr
Repeatability, fine range	sd	_	40mg+0.0006%.Rgr	_
Differential linearity deviation	sd	√(250ng⋅Rnt)	√(250ng·Rnt)	√(200ng·Rnt)
Differential eccentric load deviation	sd	0.0012%·Rnt	0.0003%·Rnt	0.0006%·Rnt
Sensitivity offset	sd	0.0008%·Rnt	0.0008%·Rnt	0.0006%·Rnt
Minimum sample weight (according to USP)		120g+0.36%.Rgr	1200g+1.8%·Rgr	120g+0.3%.Rgr
Minimum sample weight (according to USP), fine range		-	120g+1.8%·Rgr	-
Minimum sample weight (U=1%, k=2)		8g+0.024%·Rgr	80g+0.12%.Rgr	8g+0.02%·Rgr
Minimum sample weight (U=1%, k=2), fine range		_	8g+0.12%.Rgr	-
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		9.5 kg	9.5 kg	9.5 kg
Number of built-in reference weights		2	1	2
Weights for routine testing			•	•
OIML Single Weights		500 g F2	500 g F2	1000 g F2
	Small	#11118203	#11118203	#11118204
OIML Single Weights		10000 g F2	10000 g F2	20000 g F2
	Large	#11118211	#11118211	#11118212

sd = Standard deviation

Rnt = a =

Net weight (sample weight) Year (annum)

Rgr = Gross weight

¹⁾ In the temperature range 10...30 °C

		XP20000M*
Limit values		
Maximum capacity		20.1 kg
Readability		1000 mg
Tare range (fromto)		0 20.1 kg
Repeatability (at nominal load)	sd	600 mg
Linearity deviation		600 mg
Eccentricity deviation (test load)		1000 mg (10 kg)
Sensitivity offset (test weight)		800 mg (20 kg)
Sensitivity temperature drift 1)		0.001%/°C
Sensitivity stability		0.005%/a
Typical values		
Repeatability	sd	400 mg
Linearity deviation		400 mg
Eccentric deviation (test load)		600 mg (10 kg)
Sensitivity offset (test load)		240 mg (20 kg)
Minimum sample weight (according to USP)		1200 g
Minimum sample weight (U=1%, k=2)		80 g
Settling time		1 s
Dimensions		
Balance dimensions (WxDxH)		240x419x110 mm
Weighing pan dimensions		237x237 mm (WxD)
Typical uncertainties and supplementary data	1	
Repeatability	sd	400mg+0.0005%·Rgr
Differential linearity deviation	sd	√(2ug·Rnt)

		XP20000M*	
Differential eccentric load deviation	sd	0.003%·Rnt	
Sensitivity offset	sd	0.0006%·Rnt	
Minimum sample weight (according to USP)		1200g+1.5%.Rgr	
Minimum sample weight (U=1%, k=2)		80g+0.1%·Rgr	
Interface update rate		23 /s	
Weight of balance		9.5 kg	
Number of built-in reference weights		2	
Weights for routine testing			
OIML Single Weights		1000 g F2	
	Small	#11118204	
OIML Single Weights		20000 g F2	
	Large	#11118212	
sd = Standard deviation		Rnt	=
Rgr = Gross weight		a	=

¹⁾ In the temperature range 10...30 °C

* No certified models available

The precision balances with "M" platform are no longer available.

7.4.6 Balances with Readability of 0.1 g / 1 g, "L" Platform

		XP8001L	XP16001L	XP16000L
Limit values		•		-
Maximum capacity		8.1 kg	16.1 kg	16.1 kg
Readability		100 mg	100 mg	1000 mg
Tare range (fromto)		0 8.1 kg	0 16.1 kg	0 16.1 kg
Repeatability (at nominal load)	sd	80 mg	80 mg	600 mg
Linearity deviation		200 mg	200 mg	600 mg
Eccentricity deviation (test load)		300 mg (5 kg)	300 mg (5 kg)	1000 mg (5 kg)
Sensitivity offset (test weight)		640 mg (8 kg)	800 mg (16 kg)	1280 mg (16 kg)
Sensitivity temperature drift 1)		0.0015%/°C	0.0015%/°C	0.0015%/°C
Sensitivity stability		0.005%/a	0.005%/a	0.005%/a
Typical values				
Repeatability	sd	40 mg	40 mg	400 mg
Linearity deviation		130 mg	130 mg	400 mg
Eccentric deviation (test load)		200 mg (5 kg)	200 mg (5 kg)	600 mg (5 kg)
Sensitivity offset (test load)		190 mg (8 kg)	260 mg (16 kg)	380 mg (16 kg)
Minimum sample weight (according to USP)		120 g	120 g	1200 g
Minimum sample weight (U=1%, k=2)		8 g	8 g	80 g
Settling time		1.5 s	1.5 s	1.2 s
Dimensions			·	
Balance dimensions (WxDxH)		360x425x130 mm	360x425x130 mm	360x425x130 mm
Weighing pan dimensions		360x280 mm (WxD)	360x280 mm (WxD)	360x280 mm (WxD)
Typical uncertainties and supplementary data			- I	- t
Repeatability	sd	40mg+0.00025%·Rgr	40mg+0.00012%.Rgr	400mg+0.0006%·Rgr
Differential linearity deviation	sd	√(500ng⋅Rnt)	√(250ng·Rnt)	√(2.5ug·Rnt)
Differential eccentric load deviation	sd	0.002%·Rnt	0.002%·Rnt	0.006%·Rnt
Sensitivity offset	sd	0.0012%·Rnt	0.0008%·Rnt	0.0012%·Rnt
Minimum sample weight (according to USP)		120g+0.75%·Rgr	120g+0.36%.Rgr	1200g+1.8%.Rgr
Minimum sample weight (U=1%, k=2)		8g+0.05%·Rgr	8g+0.024%.Rgr	80g+0.12%·Rgr
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		12.4 kg	12.4 kg	12.4 kg
Number of built-in reference weights		2	2	2

	XP8001L	XP16001L	XP16000L			
Weights for routine testing						
OIML Single Weights	200 g F2	500 g F2	500 g F2			
Small	#11118202	#11118203	#11118203			
OIML Single Weights	5000 g F2	10000 g F2	10000 g F2			
Large	#11118206	#11118211	#11118211			
sd = Standard deviation	Rnt =	Net weight (sample weight)				

α

= Net weight (sample weight) Year (annum) =

Rgr Gross weight =

1)

In the temperature range 10...30 °C

		XP32001L	XP32001LDR	XP32000L
Limit values		1		
Maximum capacity		32.1 kg	32.1 kg	32.1 kg
Readability		100 mg	1000 mg	1 g
, Tare range (fromto)		0 32.1 kg	0 32.1 kg	0 32.1 kg
Maximum capacity, fine range		-	6.4 kg	
Readability, fine range		-	100 mg	-
Repeatability (at nominal load)	sd	80 mg	600 mg	0.6 g
Repeatability, fine range (at nominal load)	sd	-	100 mg	-
Linearity deviation		300 mg	300 mg	0.6 g
Eccentricity deviation (test load)		300 mg (10 kg)	1000 mg (10 kg)	1 g (10 kg)
Sensitivity offset (test weight)		960 mg (32 kg)	960 mg (32 kg)	1.92 g (32 kg)
Sensitivity temperature drift 1)		0.001%/°C	0.001%/°C	0.0015%/°C
Sensitivity stability		0.003%/a	0.003%/a	0.005%/a
Typical values		1	1	1
Repeatability	sd	40 mg	400 mg	0.4 g
Repeatability, fine range	sd	-	40 mg	-
Linearity deviation		200 mg	200 mg	0.4 g
Eccentric deviation (test load)		200 mg (10 kg)	600 mg (10 kg)	0.6 g (10 kg)
Sensitivity offset (test load)		320 mg (32 kg)	320 mg (32 kg)	0.65 g (32 kg)
Minimum sample weight (according to USP)		120 g	1200 g	1200 g
Minimum sample weight (according to USP), fine range		_	120 g	-
Minimum sample weight (U=1%, k=2)		8 g	80 g	80 g
Minimum sample weight (U=1%, k=2), fine range		-	8 g	-
Settling time		1.5 s	1.5 s	1.2 s
Dimensions		1	·	•
Balance dimensions (WxDxH)		360x425x130 mm	360x425x130 mm	360x425x130 mm
Weighing pan dimensions		360x280 mm (WxD)	360x280 mm (WxD)	360x280 mm (WxD)
Typical uncertainties and supplementary data			•	•
Repeatability	sd	40mg+0.00006%.Rgr	400mg+0.0003%·Rgr	0.4g+0.0003%·Rgr
Repeatability, fine range	sd	-	40mg+0.0003%.Rgr	-
Differential linearity deviation	sd	√(300ng⋅Rnt)	√(300ng·Rnt)	√(1.2ug·Rnt)
Differential eccentric load deviation	sd	0.001%·Rnt	0.003%·Rnt	0.003%·Rnt
Sensitivity offset	sd	0.0005%·Rnt	0.0005%·Rnt	0.001%·Rnt
Minimum sample weight (according to USP)		120g+0.18%.Rgr	1200g+0.9%·Rgr	1200g+0.9%.Rgr
Minimum sample weight (according to USP), fine range		-	120g+0.9%·Rgr	-
Minimum sample weight (U=1%, k=2)		8g+0.012%.Rgr	80g+0.06%.Rgr	80g+0.06%·Rgr
Minimum sample weight (U=1%, k=2), fine range		-	8g+0.06%·Rgr	-
Interface update rate		23 /s	23 /s	23 /s
Weight of balance		12.4 kg	12.4 kg	12.4 kg
5				

2

2

2

Number of built-in reference weights

Weights for routine testing

	XP32001L	XP32001LDR	XP32000L
OIML Single Weights	1000 g F2	1000 g F2	1000 g F2
Small	#11118204	#11118204	#11118204
OIML Single Weights	20000 g F2	20000 g F2	20000 g F2
Large	#11118212	#11118212	#11118212

α

sd = Standard deviation

Rgr =

Rnt = Net weight (sample weight)

= Year (annum)

¹⁾ In the temperature range 10...30 °C

Gross weight


		XP64001L	XP64000L
Limit values			
Maximum capacity		64.1 kg	64.1 kg
Readability		100 mg	1 g
Tare range (fromto)		0 64.1 kg	0 64.1 kg
Repeatability (at nominal load)	sd	100 mg	0.6 g
Linearity deviation		500 mg	0.6 g
Eccentricity deviation (test load)		500 mg (20 kg)	1 g (20 kg)
Sensitivity offset (test weight)		1280 mg (64 kg)	1.92 g (64 kg)
Sensitivity temperature drift 1)		0.001%/°C	0.0015%/°C
Sensitivity stability		0.005%/a	0.003%/a
Typical values	1		
Repeatability	sd	40 mg	0.4 g
Linearity deviation		320 mg	0.4 g
Eccentric deviation (test load)		320 mg (20 kg)	0.6 g (20 kg)
Sensitivity offset (test load)		380 mg (64 kg)	0.65 g (64 kg)
Minimum sample weight (according to USP)		120 g	1200 g
Minimum sample weight		8 g	80 g
(U=1%, k=2)		Ũ	Ū Ū
Settling time		1.8 s	1.5 s
Dimensions	•		·
Balance dimensions (WxDxH)		362x426x131 mm	362x426x131 mm
Weighing pan dimensions		362x282 mm (WxD)	362x282 mm (WxD)
Typical uncertainties and supplementary date	a		•
Repeatability	sd	40mg+0.00006%·Rgr	0.4g+0.0003%·Rgr
Differential linearity deviation	sd	√(400ng·Rnt)	√(600ng⋅Rnt)
Differential eccentric load deviation	sd	0.0008%·Rnt	0.0015%·Rnt
Sensitivity offset	sd	0.0003%·Rnt	0.0005%·Rnt
Minimum sample weight (according to USP)		120g+0.18%.Rgr	1200g+0.9%·Rgr
Minimum sample weight (U=1%, k=2)		8g+0.012%·Rgr	80g+0.06%·Rgr
Interface update rate		23 /s	23 /s
Weight of balance		14.1 kg	14.1 kg
Number of built-in reference weights		2	2
Weights for routine testing		1	- 1
OIML Single Weights		2000 g F2	2000 g F2
	Small	#11118205	#11118205
	JIIIII		
OIML Single Weights		50000 g F2	50000 g F2
	Large	#11116659	#11116659
sd = Standard deviation		Rnt =	Net weight (sample weig
Rgr = Gross weight		a =	Year (annum)

¹⁾ In the temperature range 10...30 °C


7.5 Dimensions

Dimensions in mm.


7.5.1 Balances with Readability of 0.1 mg, "S" Platform with Draft Shield

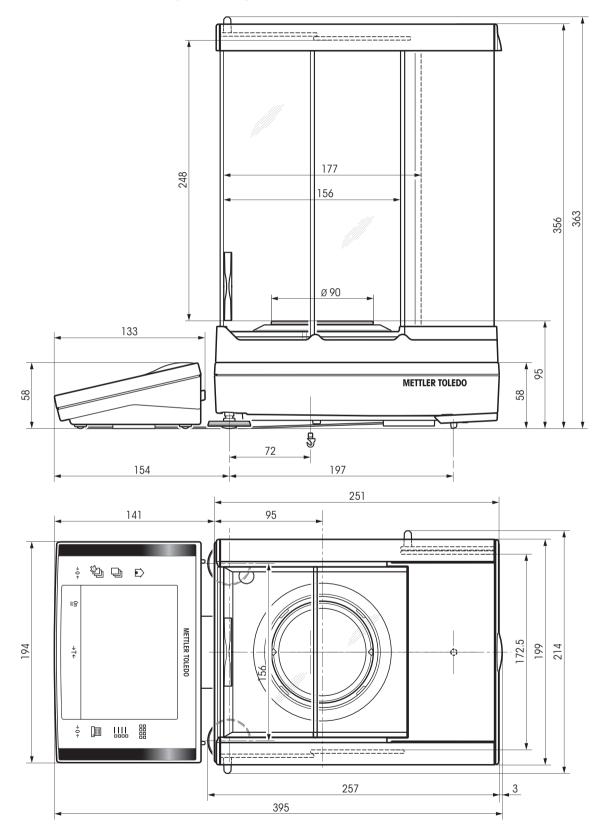


7.5.2 Balances with Readability of 1 mg, "S" Platform with Draft Shield

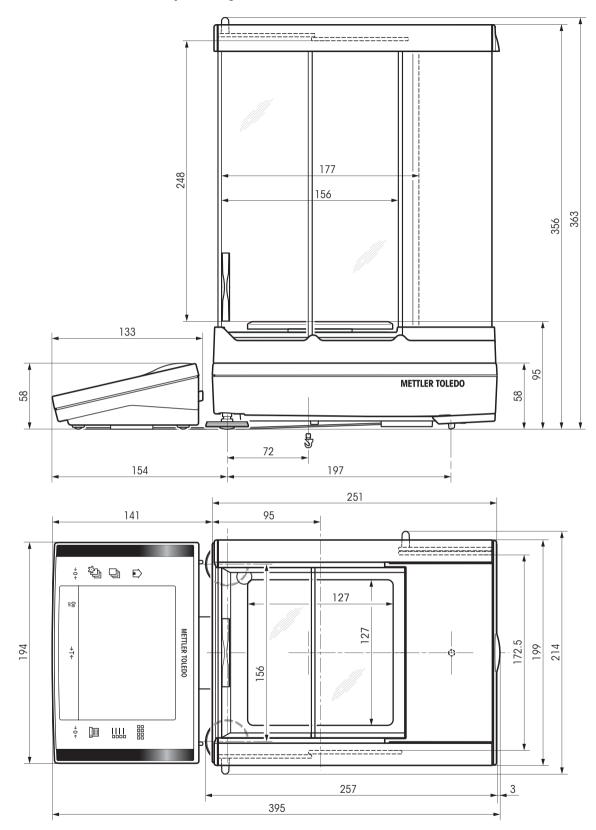


7.5.3 Balances with Readability of 10 mg, "S" Platform with Draft Shield Element

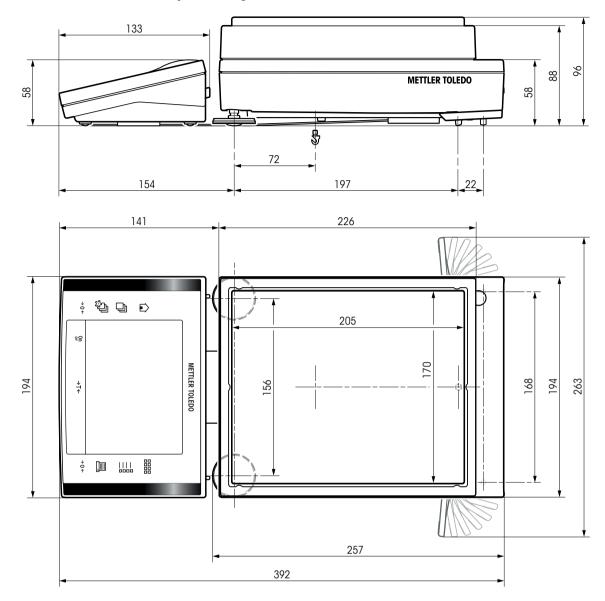


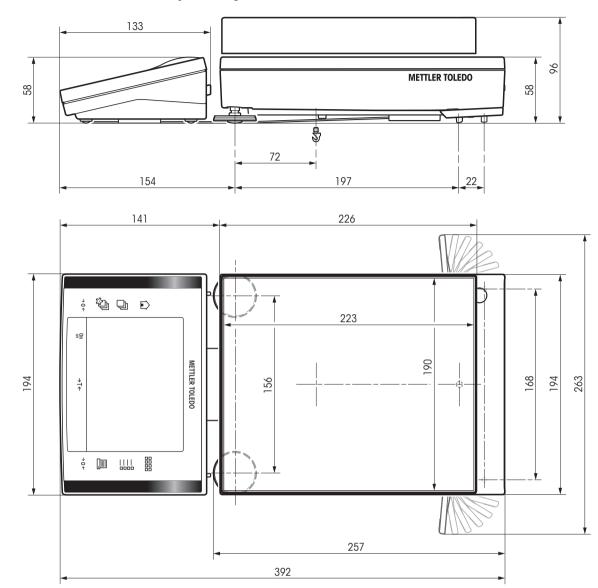


7.5.4 Balances with Readability of 0.1 g, "S" Platform


7.5.5 Balances with Readability of 10 mg / 0.1 g / 1 g, "M" Platform

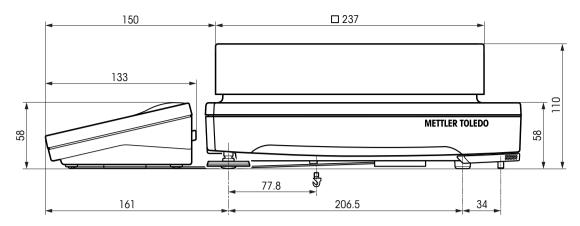

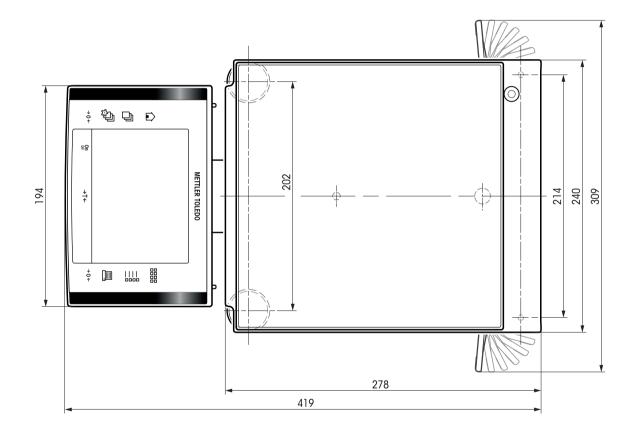

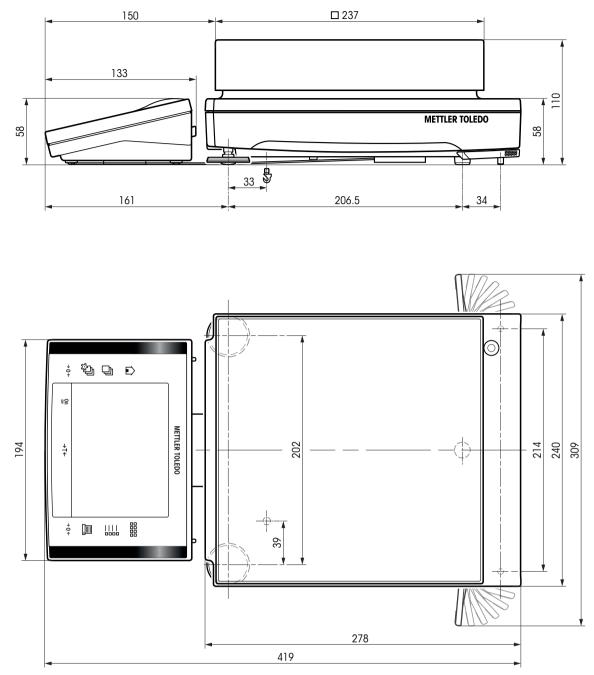

Models: XP16001M, XP16001MDR, XP20001M, XP20000M

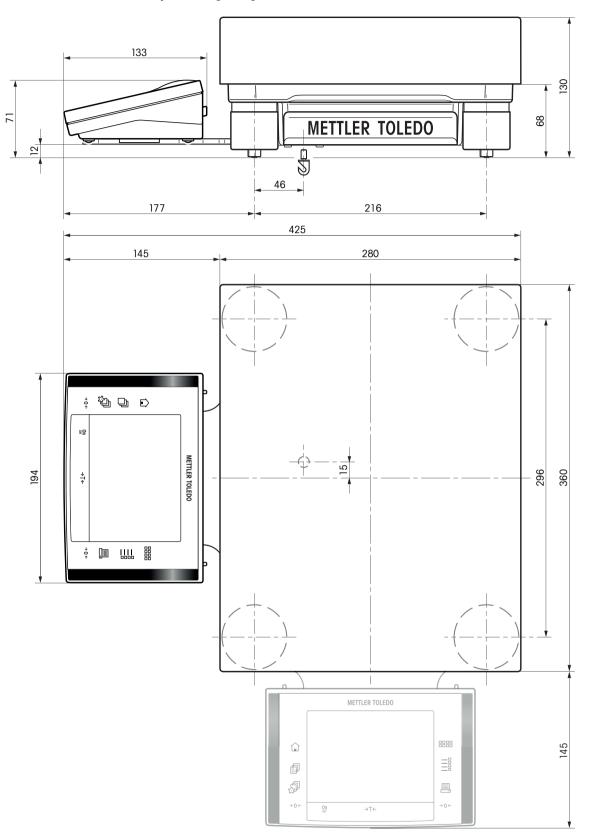

7.6 Interfaces

7.6.1 Specifications of RS232C

Interface type:	Voltage interface according to EIA RS-232C/DIN 66020 (CCITT V24/V.28)		
Max. cable length:	15 m		
Signal level:	Outputs:	Inputs:	
	+5 V +15 V (RL = 3 – 7 kΩ)	+3 V 25 V	
	–5 V –15 V (RL = 3 – 7 kΩ)	–3 V 25 V	
Connector:	Sub-D, 9-pole, female		
Operating mode:	Full duplex		
Transmission mode:	Bit-serial, asynchronous		
Transmission code:	ASCII		
Baud rates:	600, 1200, 2400, 4800, 9600, 192	00, 384001) (firmware selectable)	
Bits/parity:	7-bit/even, 7-bit/odd, 7-bit/none, 8-bit	/none (firmware selectable)	
Stop bits:	1 stop bit		
Handshake:	None, XON/XOFF, RTS/CTS (firmware s	selectable)	
End-of-line:	<cr><lf>, <cr>, <lf> (firmware se</lf></cr></lf></cr>	electable)	
GND Data	Pin 2: Balance transmit line (TxD)		
	Pin 3: Balance receive line (RxD)		
$\left \begin{array}{c} \left(\bigoplus \\ 5 \end{array} \right) \left(\bigoplus \\ 0 \end{array} \right) \left(\bigoplus \\ 1 \right) \right\rangle$	Pin 5: Ground signal (GND)		
	Pin 7: Clear to send (hardware handsl	hake) (CTS)	
Handshake	Pin 8: Request to send (hardware han	dshake) (RTS)	

7.6.2 Specifications of "Aux" Connection

You can connect the METTLER TOLEDO "ErgoSens" or an external switch to sockets "Aux 1" and "Aux 2". This allows you to start functions such as taring, zeroing, printing and others.

Connection contact

External connection Connector: 3.5 mm stereo jack connector Electrical data: Max. voltage 12 V Max. current 150 mA


8 Accessories and Spare Parts

8.1 Accessories

You can increase the functionality of your balance with accessories from the METTLER TOLEDO range. The following options are available:


	Description	Part No.
Printers		
	BT-P42 printer with Bluetooth connection to instrument	11132540
	Paper roll, set of 5 pcs	00072456
	Paper roll, self-adhesive, set of 3 pcs	11600388
	Ribbon cartridge, black, set of 2 pcs	00065975
	RS-P42 printer with RS232C connection to instrument	00229265
	Paper roll, set of 5 pcs	00072456
	Paper roll, self-adhesive, set of 3 pcs	11600388
	Ribbon cartridge, black, set of 2 pcs	00065975
	RS-P25 printer with RS232C connection to instrument	11124300
	Paper roll, set of 5 pcs	00072456
	Paper roll, self-adhesive, set of 3 pcs	11600388
	Ribbon cartridge, black, set of 2 pcs	00065975
	RS-P26 printer with RS232C connection to instrument (with date and time)	11124303
	Paper roll, set of 5 pcs	00072456
	Paper roll, self-adhesive, set of 3 pcs	11600388
	Ribbon cartridge, black, set of 2 pcs	00065975
	LC-P45 application printer with additional functions	00229119
	Paper roll, set of 5 pcs	00072456
	Paper roll, self-adhesive, set of 3 pcs	11600388
	Ribbon cartridge, black, set of 2 pcs	00065975

Optional interfaces

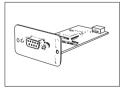

Second RS232C Interface

11132500

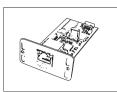

11132515



BT option: Bluetooth Interface for multipoint connection for up 11132530 to 6 Bluetooth devices



BTS option: Bluetooth Interface, single-point connection 11132535




PS/2 option: Interface for connecting commercial keyboards 11132520 and barcode readers



LocalCAN option: Interface for connection of up to 5 LC (Local-CAN) instruments



MiniMettler option: Interface MiniMettler, for connection to older	11132510
(legacy) METTLER TOLEDO systems	

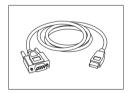

RS232 - USB converter cable – Cable with converter to connect	64088427
a balance (RS232) to a USB port	

Cables for RS232C interface


	RS9 - RS9 (m/f): connection cable for PC, length = 1 m	11101051
--	--	----------




11101052



Cables for LocalCAN Interface

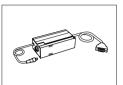

LC – RS9: Cable for connecting a PC with RS232C, 9-pin (f), 00229065 lenght = 2 m

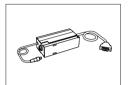

LC – RS25: Cable for connecting a printer or PC with RS232C, 00229050 25-pin (m/f), lenght = 2 m



 $LC - CL: \mbox{ Cable for connecting a device with METTLER TOLEDO } 00229130 \\ CL \mbox{ interface (5-pin), length = 2 m }$




LC - LC2: Extension cable for LocalCAN, length = 2 m	00229115
LC - LC5: Extension cable for LocalCAN, length = 5 m	00229116





LC – LCT: Cable branch (T-connector) for LocalCAN 00229118

Cables for MiniMettler Interface

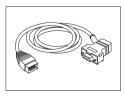

MM - RS9f: RS232C connection cable to MiniMettler interface, 00229029 length = 1.5 m

Cables for Terminal






Cable, one-sided open (2-pin)

Cable between balance and AC adapter, length = 4 m	11132037

Auxiliary displays

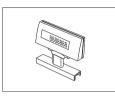

BT-BLD Bluetooth auxiliary display for table mounting, 168 mm, LCD display with backlighting	11132555
LC/RS-BLD auxiliary display on bench stand, backlit (incl. RS cable and separate AC adapter)	00224200
RS/LC-BLDS auxiliary display for table or balance mounting, 480 mm, LCD display with backlighting	11132630



RS/LC-BLDS auxiliary display for table or balance mounting,	11132630
480 mm, LCD display with backlighting	

Sensors



ErgoSens, optical sensor for hands-free operation	11132601
6 <i>i</i> 1	

LC-Switchbox



For connection of up to 3 balances with LocalCAN interface to	00229220
a printer	

Footswitches

	Footswitch with selectable function for balances (Aux 1, Aux 2)	11106741
-		





LC-FS foot switch with selectable function for balances with 00229060 LocalCAN interface

Filling-Process Control



	LV11 automatic feeder for automatic loading of small items on the balance	21900608
- A-H	Draft shield door for "Pro" Draft shield with opening for LV11	11132711
BOR	SQC14 filling process control	
	Compact instrument with printer for control of up to 16 articles	00236210
	Compact instrument with printer for control of up to 60 articles	00236211
Universal AntiStatic K	it	
11 -	Universal antistatic kit complete U-shaped, including electrode and power supply	11107767
10	Optional: Second U-electrode*	11107764
	Optional: Point-electrode*	11107765

* Power supply for optional, second U-elektrode 11107764, or for optional point-electrode 11107766 11107765

Density determination



Kit for density determination of solids and liquids (for 1 mg and 0.1 mg models)	11132680



Sinker for density of liquids in conjunction with Density Kit	00210260
Calibrated (sinker + certificate)	00210672
Recalibrated (new certificate)	00210674
Calibrated thermometer with certificate	11132685



ErgoClips



ErgoClip Sieve-S for 0.1 g and 0.01 g models, "S" platform

30039096

Ó

ErgoClip Sieve-L for "L" platform

30005760


Anti-theft devices



Steel cable

11600361

Barcode Reader

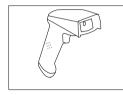

RS232C Barcode Reader		21901297
The following accessories ar	e needed for operation (not included):	
	RS232 F cable	21901305
	Null modem adapter	21900924
Plus one of the following:	AC adapter 5 V for EU	21901370
	AC adapter 5 V for US	21901372
	AC adapter 5 V for GB	21901371
	AC adapter 5 V for AU	21901370
		+ 71209966
RS232C Barcode Reader –	Cordless	21901299
The following accessories ar	e needed for operation (not included):	
	Cradle	21901300
	RS232 F cable	21901305
	Null modem adapter	21900924
Plus one of the following:	AC adapter 12 V for EU	21901373
	AC adapter 12 V for US	21901375

AC adapter 12 V for GB

AC adapter 12 V for AU



PS/2	Barcode	Reader,	without	cable
------	---------	---------	---------	-------

PS/2 wedge single cable

21901374

21901373 + 71209966

21901297

21901307



PS/2Y Barcode Reader, without cable

PS/2 wedge twin (Y) cable

21901297 21901308

Transport Cases



Transport case for precision balances for 10 mg and 0.1 g	11132595
models, "S" platform	

Protective Covers

Protective cover for XP terminal	11132570
Protective cover for weighing platform for 10 mg / 0.1 g mod- els, "S" platform	11133034
Protective cover for weighing platform, "M" platform	11132574
Protective cover for weighing pan with draft shield element for 10 mg models, "S" platform	30034950
Protective cover for weighing pan for 0.1 g models, "S" plat- form	30034949




Protective cover for weighing pan, "L" platform	30034951

Dust covers

Dust cover for 0.1 g and 0.01 g models, "S" platform

30035839

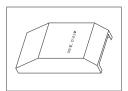

Software

LabX

•••••••		
LabX	LabX Software for One Click [™] Weighing Solutions Enables you to perform One Click [™] Standard Preparation, One Click [™] Loss on Drying, One Click [™] Sieve Analysis and many other applications. Simply start the method with the One Click [™] shortcut on the balance touchscreen. LabX guides you step-by-step through the SOP on the balance, performs your calculations automati- cally, and takes care of saving all your data. The complete solution can be tailored to match your process requirements. Visit www.mt.com/one-click-weighing for more information Freeweigh.Net	on request 21900895
M/-1-11		
Weighing pans	MPS (Magnetic Protection Shield) weighing pan for 0.1 g models 190 x 223 mm, "S" platform	11132625
	MPS (Magnetic Protection Shield) weighing pan for 10 mg models 170 x 205 mm, "S" platform	11132626
Ling V	Weighing pan 190 x 223 mm, incl. pan support, "S" platform	11132655
_/	Weighing pan 170 x 205 mm, incl. pan support and and draft shield element, "S" platform	11132660
	Dynamic weighing pan for 0.1 g and 10 mg models, with 4 litre bowl and contact mat, "S" platform	11132657

Draft shields



"MagicCube" Draft shield glass free for 0.1 mg and 1 mg mo	d- 11131650
els, usable height 175 mm, "S" platform	

11131652

"Pro" Draft shield glass free for 0.1 mg and 1 mg models (food industry), usable height 248 mm, "S" platform



Simple Draft shield glass free for 0.1 g and 10 mg models, usable height 175 mm, "S" platform	11131653
For 0.1 g model the weighing pan set must be ordered addi- tionally	11132660
Draft shield to cover the entire balance "XP-W12"	11134430



	1
300 x 450 x 450 mm (W x D x H), "S" and "M" platform	



Draft shield to cover the entire balance "XP-W64"	11134470
550 x 470 x 580 mm (W x D x H), "M" and "L" platform	

Weighing Below the Balance



Hook for the hanger (depending on the model), "M" and "L"	11132565
	11102000
platform	

Various



Terminal stand for placement of the terminal 30 cm above the weighing pan, "S" and "M" platform	11132636
Terminal stand for placement of the terminal 30 cm above the weighing pan, "L" platform	11132653







AC/DC adapter (without power cable) 100–240 VAC, 50/60 Hz, 0.3 A, 12 VDC 2.25 A	11107909
Power cable CH	00087920
Power cable EU	00087925
Power cable US	00088668
Power cable IT	00087457
Power cable DK	00087452
Power cable GB	00089405
Power cable AU	00088751
Power cable ZA	00089728
Power cable BR	30015268
Power cable JL	00225297
Power cable IN	11600569
Power cable JP	11107881
Power cable TH, PE	11107880







8.2 Spare Parts

Various

	Pos	Descriptio	n	Part No.
Weighing pans, "S" platform				
	1	0.1 mg	Weighing pan 90 mm	11133064
	2	1 mg	Weighing pan 127 x 127 mm	11131022
	3	10 mg	Weighing pan 170 x 205 mm	11131030
3	3	0.1 g	Weighing pan 190 x 223 mm	11131031
	4	10 mg	Draft shield element	11131040
		<u> </u>	, "M" platform	
4	3	all	Weighing pan 237 x 237 mm	11131173
	Weig	Weighing pans, "L" platform		
	3	≤ 32 kg	Weighing pan 280 x 360 mm	00239105
6	3	64 kg	Weighing pan 280 x 360 mm	11102124
	Varie	ous		
	5	≥ 10 mg	Pan support, "S" and "M" platform	11131029
	5		Pan support, "L" platform	00239104
	6		Foot screw, "S" and "M" platform	11106323
		Holder for	XP terminal, "S" platform	11131020
		Holder for	XP terminal, "M" platform	11131184
		XP Termin	al complete with firmware	11130690

Draft Shield, "S" platform (0.1 mg + 1 mg)

	Pos	Description	Part No.
3	1	Door front left	11133080
	2	Door back left	11133079
	3	Door top front	11133082
2	4	Door top back	11133081
	5	Door back right	11133077
	6		11133078
	7	Draft shiel element, 0.1 mg	11131531
	8	Ring seal, 0.1 mg	11131551

Packaging



Pos	Description	Part No.
"S" t	alances, Type XPxx4S	
	Packaging	11133052
	Export box	11132834
	Packaging Draft shield	11133054
	Export carton Draft shield	11132867
"S" t	alances, Type XPxx3S	
	Packaging	11133048
	Export box	11132834
	Packaging Draft shield	11133054
	Export carton Draft shield	11132867
"S" t	alances, Type XPxx2S	
	Packaging	11133046
	Export box	11132839
"S" t	alances, Type XPxx1S	
	Packaging	11133047
	Export box	11132839
"M"	balances, Type XPxxxM	
	Packaging	11133055
	Export box	11132879
"L" b	alances, Type XPxxxxL	
	Packaging	11133057
	Export box	11132912

9 Appendix

9.1 MT-SICS Interface Commands and Functions

Many of the instruments and balances used have to be capable of integration in a complex computer or data acquisition system.

To enable you to integrate balances in your system in a simple manner and utilize their capabilities to the full, most balance functions are also available as appropriate commands via the data interface.

All new METTLER TOLEDO balances launched on the market support the standardized command set "METTLER TOLEDO Standard Interface Command Set" (MT-SICS). The commands available depending on the functionality of the balance.

For further information please refer to the Reference Manual MT-SICS downloadable from the Internet under

▶<u>www.mt.com/xp-precision</u>

9.2 Procedure for Certified Balances

Preface

Certified balances are subject to the national, legal requirements of "non-automatic balances".

Switching on the balance

- Switching on
 - Immediately after being switched on, the balance displays 0.000.. g.
 - The balance is always started up with the "Factory setting" unit.

• Switch-on range

- At maximum 20% of the type load, otherwise overload is displayed (OIML R76 4.5.1).
- Stored value as switch-on zero point
 - It is not permissible to use a stored value as a switch-on zero point; the MT-SICS M35 command is not available (OIML R76 T.5.2).

Display

• Display of the weight value

- The "e" certification value is always shown in the display and is specified at the model designation plate (OIML R76 T.3.2.3 and 7.1.4).
- If the display increment is lower than the "e" certification value, this is variably displayed for the net, gross and weighed tare. (Graying of the digits or certification brackets) (OIML R76 T.2.5.4 and 3.4.1).
- In accordance with guidelines, the tested display increment (certification value) is never lower than 1 mg (OIML R76 T.3.4.2).
- At balances with d = 0.1 mg, the digits below 1 mg are displayed in gray. These digits in brackets are
 printed. In accordance with legal metrology requirements, this illustration does not affect the accuracy of the
 weighing results.

• Units of measurement

- The display and info unit are firmly set to g or mg (depending on the model).
- The following applies for the "Custom unit":
 - No certification brackets.
 - The following names are blocked, this applies to upper and lower case letters.
 - All official units (g, kg, ct etc.).
 - c, ca, car, cm, crt, cart, kt, gr, gra, gram, grm, k, kilo, to, ton.
 - All names with "o" which can be replaced by a zero (Oz, Ozt etc.).

• Identification of the weight display

- Gross, net, tare and other weight values are accordingly marked (OIML R76 4.6.5).
 - Net for net when a tare value has been used.
 - B or G for gross.
 - $-\ensuremath{\,\text{T}}$ for the weighed tare.
 - PT for the specified tare.
 - * or diff for the difference between the net or gross.

• Info field

• The info weight value is handled metrologically in the same way as the weight value in the main display.

Printout (OIML R76 4.6.11)

- If a tare value is entered manually (PreTare), the PreTare value is always printed along with the net value (PT 123.45 g).
- The printed weight values are identified in the same way as the weight value on the display. I.e. N, B or G, T, PT, diff or *, with differentiation.

Example:

Single-range balance.

Ν	123.4[5] g
PT	10.00 g \rightarrow for PreTare
G	133.4[5] g

DR balance with 100.00 g fine range.

Ν	80.4[0] g
Т	22.5[6] g \rightarrow for weighed tare
G	102.9[]g

Balance functions

Reset to zero

- The zero range is limited to a maximum of $\pm 2\%$ of the full load (OIML R76 4.5.1).
- Tare
 - No negative tare values are permitted.
 - Tare immediate (TI) is not permitted, the MT-SICS TI command is not available (OIML R76 4.6.4).
- 1/xd
 - e = d

The 1/xd switchover is not permitted (OIML R76 3.1.2).

• e = 10d

This is only permitted in the case of the 1/10d switchover.

• e = 100d

Only the 1/10d and 1/100d switchover are permitted.

10	Inde	Index				
	Α					
		AC Adapter	33			
		AC adapter	33			
		Accessories	58			
	_	Aux Connections	57			
	В		15			
		Balance cable Balance functions	15 71			
		Below-the-balance weighing	25			
	C					
		Certified balances	70			
		Cleaning	31			
		Cleaning the draft shield Conventions	31 6			
	 D	COnventions	0			
	U	Dimensions	50			
		Display	70			
		Display field remains dark	21, 22			
		Display of the weight value	70			
		Disposal	32			
	E					
		Environmental conditions ErgoSens	33, 34 57			
	F					
		Features	5			
	G					
		GLP	5			
		Good Laboratory Practice	5			
	I					
		Identification of the weight dis-	71			
		play Info field	71			
		Interface	71			
		MT-SICS	70			
		ISO 14001	5			
		ISO 9001	5			
	L					
		Level sensor	27			
		Levelcontrol Leveling Assistent	27 27			
	м	5				
		Materials	33, 34			
		MT-SICS	70			
	Р					
		Placing the terminal on the	16			
		weighing platform Power supply	20, 33, 34			
			20, 00, 04			

	Power supply voltages			20
	Printout Protection and standards		33,	71 34
R				
	Reset to zero RS232C Interface			71 57
S				
	Safety feet Safety information Scope of delivery Self test Setting the reading angle Setting up Spare Parts Switching Off Switching On		21,	28 7 13 21 22 13 68 27 27
T				
	Tare Technical data Terminal Terminal support Transport Transporting over long distances Transporting over short dis- tances Transporting the balance	15,	33, 16, 16, 23,	17 17 23 24
U				
	Units of measurement Unpacking Unpacking the draft shield			70 13 13
W				
	Weighing platform			13

GWP[®] – Good Weighing Practice™

The global weighing guideline $\mathsf{GWP}^{\circledast}$ reduces risks associated with your weighing processes and helps to

- choose the appropriate balance
- reduce costs by optimizing testing procedures
- comply with the most common regulatory requirements

www.mt.com/GWP

www.mt.com/excellence

For more information

Mettler-Toledo AG, Laboratory & Weighing Technologies CH-8606 Greifensee, Switzerland Tel. +41 (0)44 944 22 11 Fax +41 (0)44 944 30 60 www.mt.com

Subject to technical changes. © Mettler-Toledo AG 05/2012 11781055B en

