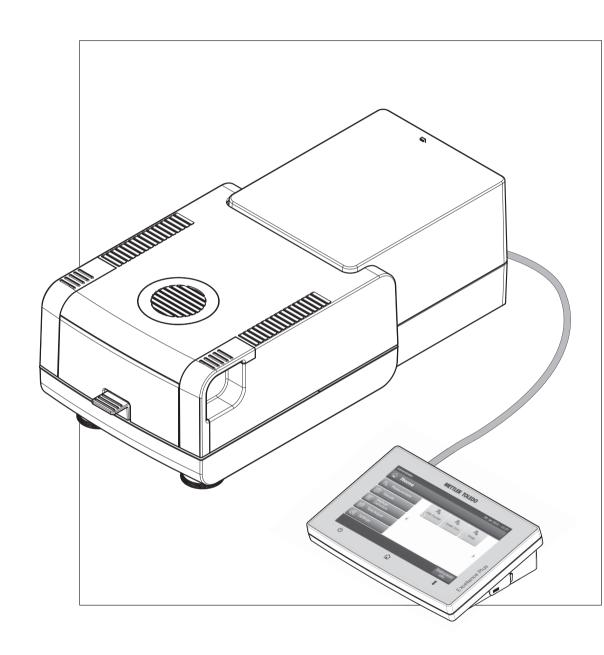
Moisture Analyzer

Excellence Plus HX204



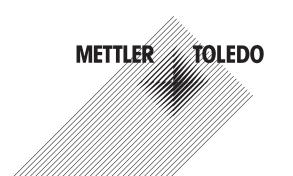


Table of contents

1	Introductio	ns		7
		1.1	Conventions and Symbols Used in these Operating Instructions	7
2	Safety Info	rmation		8
		2.1	Definition of Signal Warnings and Symbols	8
		2.2	Product Specific Safety Notes	8
3	Design and	function		12
		3.1	Overview	12
		3.1.1	Drying Unit	12
		3.1.2	Terminal	13
		3.1.3	Keys on the Terminal	15
		3.2 3.2.1	User Interface User "Home" Screen	16 17
		3.2.2	Basic Elements on Touch Screen	18
		3.2.3	Icon and Buttons	18
		3.2.3.1	Status Message Icons	18
		3.2.3.2	Title Bar Icons	18
		3.2.3.3 3.2.4	Action Bar Buttons Input Dialogs	18 19
		3.2.4.1	Entering Text and Numbers	19
		3.2.4.2	Entering Numerical Values	19
		3.2.4.3	Changing Date and Time	20
		3.2.5	Lists and Tables	21
		3.2.6 3.2.7	Work Screen Graphics View	21 22
4	Installation		·	23
		4.1	Scope of Delivery	23
		4.2	Location	23
		4.3	Opening and Closing Sample Chamber	24
		4.4	Connecting Instrument	24
		4.5	Setting up Drying Unit	25
		4.6	Leveling Drying Unit	26
		4.7	Anti-theft Device	27
		4.7	Setting Reading Angle of Terminal	27
		4.9	Setting Date and Time	27
		4.10	Adjustment	28
5	My First M		·	29
	<u> </u>	Gusureniei	"	
6	Settings	0.1	H . D .	32
		6.1 6.1.1	User Preferences	33 33
		6.1.1	Languages Screen	33
		6.1.3	Password	33
		6.2	System Settings	34
		6.2.1	Regional Settings	34
		6.2.2	Peripherals	34

		6.2.3 6.2.3.1 6.2.3.2 6.2.4	Default Settings Default Languages Default Screen Settings Touch Screen Adjustment	35 35 36 36
		6.3 6.3.1 6.3.2	Applications Settings Identifications Management Print and Output Management	38 38 38
		6.4 6.4.1.1 6.4.1.2 6.4.1.3 6.4.1.4 6.4.1.5 6.5 6.5.1 6.5.2		40 40 40 41 42 42 43 44 44 44
		6.5.3 6.6 6.6.1 6.6.2 6.6.3 6.6.4	Account Policies System- and Data Management Export / Import Back up / Restore Reset Update	47 48 48 49 50 50
7	Test/Adjust	7.1 7.1.1 7.1.2	Adjustments Weight Adjustment - Internal Weight Adjustment - External	52 52 53 53
		7.1.3 7.1.4 7.2 7.2.1 7.2.2 7.2.3 7.2.4 7.3	Temperature Adjustment Adjustment Printout Examples Tests Weight test - internal Weight test - external Temperature test Test Printout Examples History	54 56 57 57 57 57 58 59
8	Method Def	inition	·	61
		8.1 8.1.1.1 8.1.1.2 8.1.1.3 8.1.1.4 8.1.1.5 8.1.2 8.1.3 8.2 8.2.1 8.2.2	Main Measuring Parameters Settings of Drying Program Settings of Standard drying Settings of Rapid drying Settings of Gentle drying Settings of Step drying Settings of Switch-off Criterion Settings of Display Mode Settings of Start Weight Result and Values Handling Control limits Resolution	63 63 63 64 65 66 67 69 71 72 72

		8.3 8.3.1	Workflow Handling Starting Mode	75 75
		8.3.2	Preheating	75
		8.4 8.4.1	General Method Properties Method name	77 77
		8.5	Test a Method	78
9	Mogourom		resi a Mellioa	
9	Measurem		B (' ' N	80
		9.1	Performing a Measurement	80
		9.2	Work with Weighing-in Aid	82
		9.3	Using Shortcuts	82
10	Results			84
		10.1	Graphical Evaluations of Measurement Results	84
		10.2	Information on Printouts	86
		10.3	Exporting Results	87
11	Maintenan	ce		88
		11.1	Cleaning	88
		11.1.1	Sample Chamber	89
		11.1.2	Heating Module	89
		11.1.3	Fan Grill	91
		11.2	Dust Filter	91
		11.3	Replacing Power Line Fuse	92
		11.4	Disposal	92
12	Troublesho	•		93
		12.1	Error Messages	93
		12.2	Status Messages	93
		12.3	What to Do if	95
13	Technical I	Data		97
		13.1	General Technical Data	97
		13.2	Dimensions	99
14	Accessorie	s and Spa	re Parts	100
15	Appendix			104
		15.1	How to Obtain Best Results	104
		15.1.1	Measurement Principle of Halogen Moisture Analyzer	104
		15.1.2	Notes on Adjustment of Balance and Heating Module	104
		15.1.3	Optimum Sample Preparation	105
		15.1.4	Further Information on Moisture Determination	105
		15.2	Recommended Printer Settings	105
		15.3	Third Party License/Notice	106
16	Indov			107

1 Introductions

Thank you for purchasing a Halogen Moisture Analyzer from METTLER TOLEDO.

Your Moisture Analyzer is fast and reliable. It offers a high level of operating convenience and useful functions to facilitate determination of the moisture content of your samples.

Behind your instrument stands METTLER TOLEDO, a leading manufacturer of not only balances and scales for the lab and production, but also analytical measuring instruments. A customer service network covering the entire globe with well trained personnel is your service at all times, whether you are choosing accessories or require guidance for a specific application to ensure optimum utilization of your instrument.

Your Halogen Moisture Analyzer is used for determining the moisture content of almost any substance. The instrument works on the thermogravimetric principle. At the start of the measurement the Moisture Analyzer determines the weight of the sample, the sample is then quickly heated by the integral halogen heating module and the moisture vaporizes. During the drying process the instrument continually measures the weight of the sample and displays the reduction in moisture. Once drying has been completed, the moisture or solids content of your sample is displayed as the final result.

Of decisive importance in practice is the rate of heating and even heating of the surface of the sample. In comparison with conventional infrared heating or the drying oven method, for example, the halogen heating module of your instrument needs a shorter time to reach its maximum heating power. It also allows use of high temperatures, an additional factor in shortening the drying time. Uniform heating of the sample material ensures good repeatability of the drying results and makes it possible to use a smaller amount of sample.

The Moisture Analyzer conforms with all common standards and directives. It supports stipulations, work techniques and result records as demanded by all international quality assurance systems, e.g. GLP (Good Laboratory Practice), GMP (Good Manufacturing Practice). The instrument has a CE declaration of conformity and METTLER TOLEDO as the manufacturer has been awarded ISO 9001 and ISO 14001 certification. This provides you with the assurance that your capital investment is protected in the long term by a high product quality and a comprehensive service package (repairs, maintenance, servicing, adjustment service).

Finding More Information

▶ www.mt.com/moisture

Software version

These operating instructions are based on the initially installed firmware (software) version V1.0.

1.1 Conventions and Symbols Used in these Operating Instructions

Key and button designations are indicated by a picture or text in square brackets (e.g. [Save]

These symbols indicate an instruction:

- prerequisites
- ı steps
- 2 ...

2 Safety Information

2.1 Definition of Signal Warnings and Symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

Signal Words

WARNING for a hazardous situation with medium risk, possibly resulting in severe

injuries or death if not avoided.

CAUTION for a hazardous situation with low risk, resulting in damaged to the device

or the property or in losing of data or minor or medium injuries if not avoid-

ed.

Attention (no symbol)

for important information about the product.

Note (no symbol)

for useful information about the product.

Warning Symbols



General hazard



Electrical shock



Hot surface



Inflammable or explosive substance



Toxic substance



Acid / Corrosion

2.2 Product Specific Safety Notes

General Safety Information

Your instrument meets the state of the art technology and complies with all recognized safety rules, however, certain hazards may arise in extraneous cirmumstances. Do not open the housing of the instrument: It does not contain any parts which can be maintained, repaired or replaced by the user. If you ever have problems with your instrument, contact your authorized METTLER TOLEDO dealer or service representative.

Always operate and use your instrument only in accordance with the instructions contained in this manual. The instructions for setting up your new instrument must be strictly observed.

If the instrument is not used according to these Operating Instructions, protection of the instrument may be impaired and METTLER TOLEDO assumes no liability.

Intended Use

Your Moisture Analyzer is used for determining the moisture in samples. Use the instrument exclusively for this purposes. Any other type of use and operation beyond the limits of technical specifications without written consent from Mettler-Toledo AG, is considered as not intended.

Moisture determination applications must be optimized and validated by the user according to local regulations. Application-specific data provided by METTLER TOLEDO is intended for guidance only.



It is not permitted to use the instrument in explosive atmosphere of gases, steam, fog, dust and flammable dust (hazardous environments).

Staff Safety

The moisture Analyzer may be operated only by trained personnel who are familiar with the properties of the samples used and with the handling of the instrument.

In order to use the instrument, you must have read and understood the operating instructions. Keep the operating instructions for further reference.

Never make any modifications or constructional alterations to the instrument and use only original spare parts and optional equipment from METTLER TOLEDO.

Protective Clothing

It is advisable to wear protective clothing in the laboratory when working with the instrument.



A lab coat should be worn.



A suitable eye protection such as goggles should be worn.



Use appropriate gloves when handling chemicals or hazardous substances, checking their integrity before use.

Safety Notes



WARNING

Risk of electric shock

Your instrument is supplied with a 3-pin power cable with an equipment grounding conductor. Only extension cables which meet this relevant standards and also have an equipment grounding conductor may be used. Intentional disconnection of the equipment grounding conductor is prohibited.

CAUTION



The Halogen Moisture Analyzer works with heat!

- a) Ensure sufficient free space around the instrument to avoid heat accumulation and overheating (approx. 1 m free space above the heating module).
- b) The vent over the sample must never be covered, plugged, taped over or tampered with in any other way.
- c) Do not place any combustible materials on, under or next to the instrument since the area around the heating module may be hot.
- d) Exercise caution when removing the sample. The sample itself, the sample chamber, the draft shield and any sample vessels used may still be very hot.
- e) During operation, you should never open the heating module itself as the ring-shaped heating reflector or its protective glass can reach 400 °C! If you have to open the heating module e.g. for maintenance, disconnect the instrument from the power supply and wait until the heating module has cooled down completely.
- f) No modifications must be made within the heating module. It is particularly dangerous to bend any components or remove them or to make any other changes.

Certain samples require special care!

With certain types of samples, there is a possibility of danger to personnel or damage of property. Please note that the user always has the responsibility and liability for damage caused by use of any types of samples!



CAUTION

Fire or Explosion

- Flammable or explosive substances.
- Substances containing solvents.
- Substances which evolve flammable or explosive gases or vapors when heated.
- a) In cases of doubt, perform a careful risk analysis.
- b) Work at a drying temperature that is low enough to prevent the formation of flames or an explosion.
- c) Wear protective goggles.
- d) Work with small amounts of sample.
- e) Never leave the instrument unattended!



WARNING

Substances which contain toxic or caustic components

Toxic gases produced during drying could cause irritations (eyes, Skin, breathing), illness or death.

Such substances may be dried only in a fume cupboard.

CAUTION



Corrosion

Substances which evolve corrosive vapors when heated (e.g. acids).

 Work with small amounts of samples as the vapor can condense on cooler housing parts and cause corrosion.


3 Design and Function

In order to work with your Moisture Analyzer, you need to familiarize yourself with the instrument layout, drying unit, graphical user interface of the touch screen as well as the terminal.

3.1 Overview

3.1.1 Drying Unit

Description of the components and the external connectors on the back of the drying unit (Weighing unit with heating module).

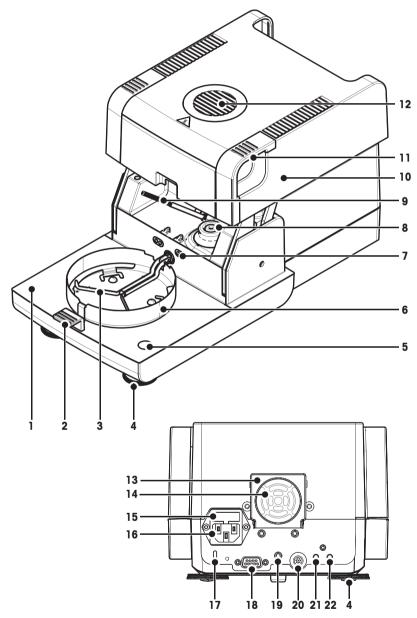


- 1 Sample chamber
- 2 Sample pan handler
- 3 Sample pan holder
- 4 Leveling screw
- **5** Level indicator / Level sensor
- 6 Draft shield element
- 7 Contacts for temperature adjustment kit(optional)
- 8 Power contacts for halogen lamp
- 9 Temperature sensor
- **10** Heating module (with thermal overload protection)
- 11 Handles for emergency opening the sample chamber
- 12 Inspection window and vent
- **13** Fan
- 14 Dust filter (optional)
- 15 Power line fuse
- **16** Power supply socket
- 17 Kensington slot for antitheft purposes
- **18** Humidity sensor connection (n.a.)
- **19** Fastening for terminal stand (optional)
- **20** System connection (terminal cable)
- 21 Aux 2 connection (n.a.)
- 22 Aux 1 connection (n.a.)

Connectivity

System connection	Interface for the communication between terminal and drying unit.
Power supply socket	Power supply inlet socket for the country-specific power cable.
	The socket has a built in fuse compartment containing a replacement fuse.

3.1.2 Terminal

Description of the connections at the terminal (Input and output unit with interfaces).

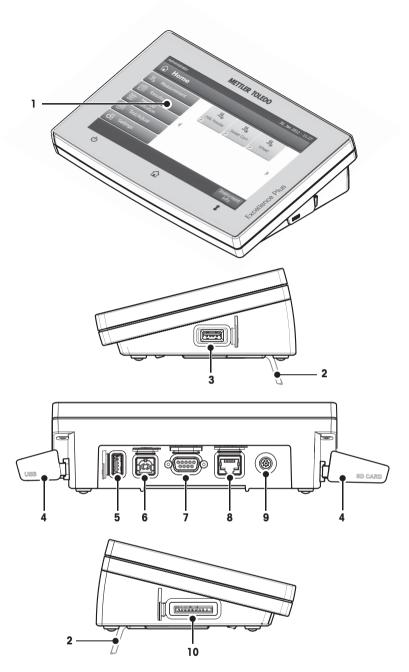


- 1 WVGA color display (touch screen)
- 2 Foot
- 3 USB Host connection
- 4 Faceplate
- 5 USB Host connection
- **6** USB Device connection (n.a.)
- 7 RS232C connection
- **8** Ethernet RJ45 connection (n.a.)
- **9** System connection (terminal cable)
- **10** Memory Card slot for SD Card

Connectivity

••		
System connection	Interface for the communication between terminal and drying unit.	
RS232C connection	RS232C interface for connecting the instrument to a peripheral device (e.g. printer or a PC with a 9-pin male connector).	
USB Host connection	USB interface for connecting a peripheral device to the instrument (e.g. Memory Stick). Type A Socket.	
Memory Card slot	The memory card slot fits Secure Digital Memory Cards. SD and SDHC types as well as sizes up to 32 GB are supported. You can use it as an external storage device e.g. to store measured results.	
	Note SDXC types are not supported.	

3.1.3 Keys on the Terminal

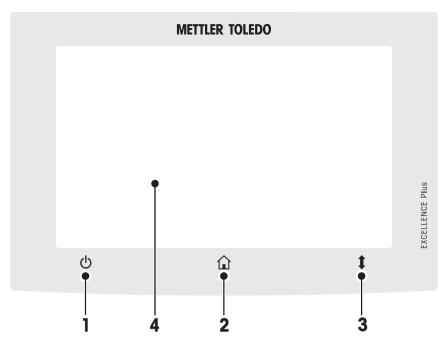

		Name	Explanation
1	داء	ON/OFF	To switch on or off (standby operation) the instrument.
			Note Do not disconnect the instrument from the power supply except if you will not be using the instrument for an extended period.
2		Home	To return from any menu level, or other window to the user home screen.
3	‡	Open/Close	To open or close the sample chamber motorized. Attention Do not open the sample chamber manually. Only open manually in emergency situations. In all other cases use always the [‡] key. See Opening and Closing Sample Chamber (page 24)
4		Touch screen	

3.2 User Interface

The colored touch screen is a touch-sensitive WVGA monitor. The touch screen not only displays information, it also allows you to enter commands by touching certain areas on its surface: You can choose the information displayed on the screen, change terminal settings or perform certain operations on the instrument.

Only those elements which are available for the current dialog appear on the display.

Buttons on the Touch Screen

Buttons are software elements on the touch screen (Soft keys).




CAUTION

Do not touch the touch screen with pointed or sharp objects!

This may damage the touch screen.

3.2.1 User "Home" Screen

The user **Home** screen is the main screen and appears after startup or login of the instrument. All screens of the user interface can be accessed from the home screen. Returning to the **Home** screen is possible from all other screens of the user interface by pressing the $[\widehat{\ }]$ key or tapping the $[\mathbf{Home}]$ button.

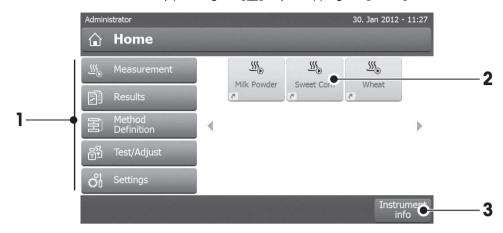


	Name	Explanation
1	Measurement	Start measurement
		Prerequisite: Method is defined and all necessary settings are made.
		For further information see My First Measurement (page 29) and Measurement (page 80).
	Results	Show results
		Under this menu item, all measurement results can be displayed, printed or exported. For further information see Results (page 84).
	Method Definition	Define, edit or delete a method
		All actions using in a drying method, can be defined under this menu item. Description of the menu item see Method Definition (page 61).
	Test/Adjust	Adjust or test the instrument
		Under this menu item, the integrated balance and the heating module can be adjusted or tested.
		Description of the menu item see Test/Adjust (page 52).
	Settings	Define general settings
		All settings of the system can be defined, e.g, system settings, user management or user preferences. These settings are usually made during installation of the instrument. Description of the menu item see Settings (page 32)
2		User shortcuts
_		User-specific shortcuts for frequently used methods. Shortcuts are saved in the
		user profile. How to create shortcuts see Using Shortcuts (page 82).
3	Instrument info	Shows general info about the instrument like type, serial number, software version.

3.2.2 Basic Elements on Touch Screen

The graphical user interface consists of the following basic elements.

Screen Areas and their Elements

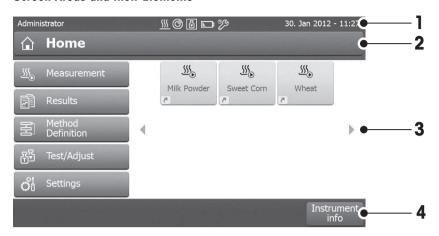

	Name	Explanation
1	Status Bar	The status bar contains the user name, status icons as well as date and time. See Status Message Icons (page 18).
2	Title Bar	The title bar contains elements for the user's orientation and information, as well as for special functions. See Title Bar Icons (page 18).
3	Content Area	The content area is the main working area of the menus and the applications, content depends on application or performed action. The content area also shows graphic illustrations such as the curve diagram of a data set (e.g. drying curve).
4	Action Bar	The action bar contains action buttons for triggering the actions which are required and available in the current dialog (e.g. Back , -> 0/T<- , Print , Save , Delete , OK). See Action Bar Buttons (page 18).

3.2.3 Icon and Buttons

3.2.3.1 Status Message Icons

Status messages are displayed by means of small icons in the status bar. For the most important messages **see** Status Messages (page 93).

By tapping the symbols, the explanation of the symbols will appear in a separate window.

3.2.3.2 Title Bar Icons

The title bar icons only appear when they are active. The icons which appear in the title bar, are listed with their function below:

	Explanation
[5]	Create a shortcut and add it to your user home screen. Edit or remove an existing shortcut
	see Using Shortcuts (page 82).

3.2.3.3 Action Bar Buttons

The action bar contains action buttons for triggering any of the actions which are required and available in the current dialog (e.g. **Back**, **->0/T<-**, **Print**, **Save**, **Delete**, **OK**). There are up to six action buttons available for a work step. The functions of these buttons are self explanatory and therefore not discussed further.

3.2.4 Input Dialogs

3.2.4.1 Entering Text and Numbers

The keyboard dialog allows you to enter characters including letters, numbers and a variety of special characters. The country-specific layout of the keyboard is in accordance with the selected language, **see** Regional Settings (page 34).



	Name	Explanation
1	Input field	
2	Explanation field	The maximum number of characters allowed is indicated.
3	Shiff	You can enter lower and upper case letters by toggling 😭. Switch the keyboard
		by tapping the corresponding button in the action bar, for entering numbers, spe-
		cial characters or letters

Confirm with [OK] (To exit keyboard tap [Cancel]).

	Function
핲	Shift-Lock
\boxtimes	Delete last character
(Cursor left
\rightarrow	Cursor right

Note

It is also possible to place cursor directly in input field by tapping the respective position.

3.2.4.2 Entering Numerical Values

The numerical keypad dialog allows you to enter numerical values.



	Name	Explanation
1	Switch	If a switch appears, located on the left side in the title bar; it must be switched on
		for the entry of a value: $O = OFF$, $I = ON$
2	Input field	
3	Explanation field	The valid input range is indicated.


Confirm with [OK].

To exit keypad, tap [Cancel].

	Function	
\boxtimes	Delete last character	
←	Cursor left	
\rightarrow	Cursor right	

3.2.4.3 Changing Date and Time

This dialog (Picker view) allows you to set the date and time by picking from the built-in calendar/clock.

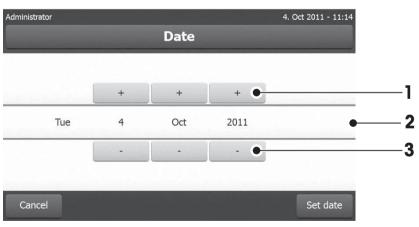

	Name	Explanation
1	Pick button	Pick button upwards (input control)
2	Explanation field	Current value
3	Pick button	Pick button downwards (input control)

- Confirm with [Set date]. (To exit without changing, tap [Cancel].)

3.2.5 Lists and Tables

Simple list view

Basic elements of a simple list including a content title. If necessary, the content area can be scrolled vertically and can be arranged in a different order.



	Name	Explanation
1	Content title	
2	Content area	
3	Arrow icon	[↓] Showing the list in ascending order.[↑] Showing the list in descending order.
		[1] Showing the list in descending order.
4	Scrollbar	Scrolling up
5		Scroller
6		Scrolling down

Further types of lists:




 $[\wedge]$

Accordion panels can be opened or closed by tapping Closable panels can be opened or closed by tapping [>].

3.2.6 Work Screen

The work screen is used for performing main tasks e.g. measurements.

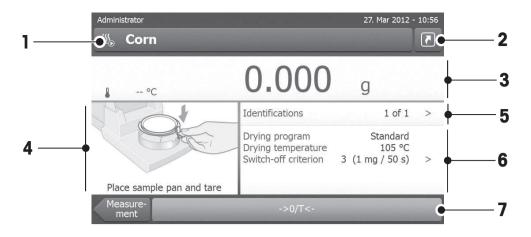

	Name	Explanation	
1	Name field	Name of current Method	
2	Shortcut button	Adding/editing a shortcut to the home screen for this method	
3	Value panel	Displays the current measured values of the work process	
4	Graphics panel	E.g. graphical illustrations of the drying curves, instructions for the user for per-	
		forming tasks, and weighing-in aid	
5	ID panel	Identifications (IDs) appear after tapping the ID panel for entering or editing val-	
		ues (comments). ID panel appears only if identification input is activated in the	
		menu.	
6	Parameter panel	Displays the parameters of the current work process. A detailed overview of the	
		method parameters appears after tapping the parameter panel.	
7	Action buttons	Depending on the current context	



Measurement in progress

3.2.7 Graphics View





Coffee Powder

5.53 %MC

Graphical view of single measurement

For further information see Graphical Evaluations of Measurement Results (page 84)

4 Installation and Commissioning

In this section you will find information how to put your new instrument into operation.

4.1 Scope of Delivery

Open the package and remove drying unit, terminal and the accessories. Check the completeness of the delivery. The following accessories are part of the standard equipment of your new Moisture Analyzer:

- 80 aluminum sample pans
- 1 sample handler
- 1 sample pan holder
- 1 specimen sample (circular, absorbent glass fibre filter)
- 1 draft shield
- 1 power cable (country specific)
- 1 Terminal cable (Interface Terminal drying unit)
- 1 Operating instructions
- 1 Application brochure «Guide to Moisture Analysis»
- 1 CE declaration of conformity

Remove the packaging from the instrument. Check the instrument for transport damage. Immediately inform your METTLER TOLEDO representative if you have any complaints or parts are missing.

Note

Store all parts of the packaging. This packaging guarantees the best possible protection for the transport of your instrument.

4.2 Location

Your Moisture Analyzer is a precision instrument. An optimum location guarantees accuracy and dependability. Make sure that the following environmental conditions are met:

- Operate the instrument only indoors and at an altitude of less than 4000 m above sea level.
- Before switching on the instrument, allow all its parts to reach room temperature (+5 to 30°C).
 Make sure that the relative humidity is between 20% and 80% and non-condensing conditions are met.
- The power plug must be easily accessible.
- Firm, horizontal location as free from vibrations as possible.
- Avoid direct sunlight.
- No excessive temperature fluctuations.
- No powerful drafts.
- Surroundings as free from dust as possible.
- Sufficient clearance around the instrument to allow warm air to dissipate.
- Sufficient distance from heat-sensitive materials in the vicinity of the instrument.









WARNING



Substances which contain toxic or caustic components

Toxic gases produced during drying could cause irritations (eyes, Skin, breathing), illness or death.

Such substances may be dried only in a fume cupboard.

4.3 Opening and Closing Sample Chamber

Open and close the sample chamber exclusively by pressing the [1] key.

Attention

Do not open the sample chamber manually.

Only manually opening the sample chamber in emergency situation:

- 1 Hold the handles on the heating module **with both hands** and pull the module firmly upwards to open the sample chamber.
- 2 Pull out the power plug.

Note

It is necessary to reboot the system, after an emergency opening.

4.4 Connecting Instrument



WARNING

Risk of electric shock

- a) Use only the 3-pin power cable with equipment grounding conductor which was supplied with your instrument.
- b) Only an 3-pin grounded outlet for connecting your instrument must be used.
- c) Only extension cables which meet this relevant standards and also have an equipment grounding conductor may be used.
- d) Intentional disconnection of the equipment grounding conductor is prohibited.

Attention

First, check whether the voltage printed on the type plate of the drying unit matches your local line voltage. If
this is not the case, on no account connect the drying unit to the power, but contact your METTLER TOLEDO
sales representative or dealer.

Two different versions of drying units with country-specific power cable are available (110 V AC or 230 V AC).

- The power plug must be always accessible.
- Before operating, check all cables for damage.
- Guide the cables so that they cannot become damaged or interfere with the measuring process.

Connecting the Terminal to the Drying Unit


- Drying unit and terminal are at the final location.
- 1 Connect one end of the supplied terminal cable to the system connection socket on the terminal, **see** Terminal (page 13).
- 2 Screw the connector firmly.
- 3 Connect the other end to the system connection socket on the drying unit, see Drying Unit (page 12).

4 Screw the connector firmly.

Connecting the Drying Unit to the Power Supply

- ► Terminal drying unit interface is connected.
- 1 Connect one end of the power cable to the power supply socket on the drying unit, **see** Drying Unit (page 12).
- 2 Connect the other end to the power line outlet.

4.5 Setting up Drying Unit

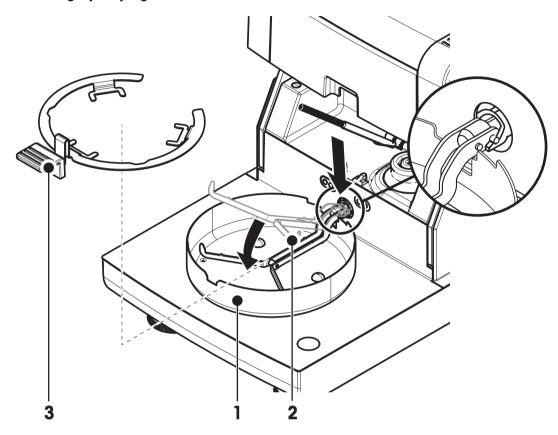

- ▶ Terminal and drying unit are connected.
- 1 Switch on the Instrument by pressing [4].
- 2 Press [1].
 - ⇒ the motorized automatic sample chamber opens.
- 3 Place the draft shield element (1). Only one position is possible.
- 4 Insert the sample pan holder (2) carefully. Check that the sample pan holder is hooked correctly.
- 5 Insert the sample pan handler (3).

Attention

The drying unit is specially designed for motorized opening or closing of the sample chamber. Do not open the sample chamber manually. Only open manually in emergency situations. In all other cases use always the [‡] key, see Keys on the Terminal (page 15).

Note

When putting into operation for the first time or if the instrument is disconnected from the power over a longer period, leave the instrument connected to the power supply for at least 5 hours to allow the built-in rechargeable battery to charge up! This battery ensures that the date and time are not lost when the instrument is dis-

connected from the power. This battery can not be replaced by the user. Please contact a METTLER TOLEDO representative.


4.6 Leveling Drying Unit


Exact horizontal positioning and stable installation are prerequisites for repeatable and accurate results. To compensate for small irregularities or inclinations (±2 %) at the location, the instrument must be leveled.

For exact horizontal positioning, the drying unit has a level indicator (level) and 2 leveling screws. When the air bubble in the level indicator is exactly in the center, the instrument is standing perfectly horizontally. To level it, proceed as follows:


- 1 Position your Moisture Analyzer at the selected location.
- 2 Turn the two leveling screws until the air bubble is in the center of the level indicator.

L = left foot R = right foot

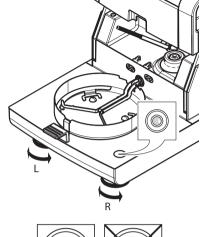

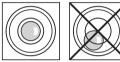

Air bubble at	"12 o'clock"	turn both feet clockwise
Air bubble at	"3 o'clock"	turn left foot clockwise, right foot counterclock- wise
Air bubble at	"6 o'clock"	turn both feet counter- clockwise
Air bubble at	"9 o'clock"	turn left foot counterclock- wise, right foot clockwise

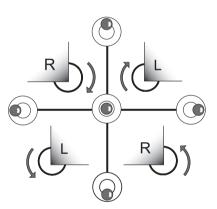


Note

The drying unit should be re-leveled each time its location is changed.

Leveling the drying unit with optional stand for terminal mounted.

Stand for terminal **see** Accessories and Spare Parts (page 100).

- The drying unit is positioned at the selected location
- 1 To loose the clamps for the safety feet of the stand, turn them outwards.
- 2 Level the drying unit as previously described.
- 3 To lock the clamps for the safety feet, turn them inwards as far as they will go.

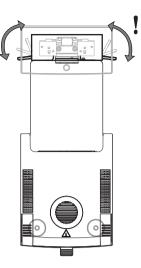

Level sensor


The built-in level sensor constantly monitors the correct horizontal alignment of the drying unit. When the level is incorrect, the status icon @ appears in the display, **see** Status Message Icons (page 18).

Note

The level sensor cannot be switched off.

4.7 Anti-theft Device

For protection against theft, your drying unit is fitted with a Kensington security slot for an anti-theft device. To find a suitable anti-theft cable **see** Accessories and Spare Parts (page 100).

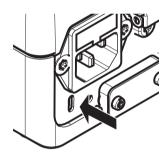

4.8 Setting Reading Angle of Terminal

- To change the reading angle, fold out the two tilt feet.



4.9 Setting Date and Time

Navigation: Home > Settings > System settings > Regional Settings

When you put your new instrument into operation for the first time, you should enter the current date and time. These settings are retained even if you disconnect your instrument from the power supply.

Note

Various languages are available for the dialog with your instrument and can be selected in the menu: **Navigation: Settings** > **User preferences**, **see** User Preferences (page 33).

Set the current date

- ▶ Regional Settings is selected.
- 1 Tap Date.
- 2 Set day, month and year.
- 3 Confirm with [Set date].

Set the current time

- ▶ Regional Settings is selected.
- 1 Tap **Time**.
- 2 Set Hours and minutes.
- 3 Confirm with [Set time].

Note

For changing the format for date and time, see Regional Settings (page 34).

4.10 Adjustment

To obtain accurate measuring results, it is necessary to adjust the integrated balance as well as the heating module.

Adjusting is necessary:

- before the instrument is used for the first time.
- at regular intervals.
- after a change of location.

The following adjustment options are possible:

- Fully automatic balance adjustment FACT.
- Balance adjustment with internal weight.
- Balance adjustment with external weight (accessory).
- Temperature adjustment with temperature adjustment kit (accessory).

For more information to perform these operations see Test/Adjust (page 52).

5 My First Measurement

After you have successfully put your new Moisture Analyzer into operation for the first time, you can immediately perform your first measurement. In doing so, you will become familiar with the instrument.

Use the supplied specimen sample (absorbent glass fiber filter) for your first measurement to determine the moisture content. During your first measurement the instrument operates with the factory settings.

Switch on

- ▶ The instrument is connected to the mains.
- To switch on, press [(1)].



Create a method

- 1 Tap Method Definition.
 - Method Definition menu appears.
- 2 To define a new method, tap [New...].
 - ⇒ The keypad appears.
- 3 Enter a name for your first method e.g. My first method.
- 4 Confirm with [**OK**].
- 5 To store your new method with factory settings, tap [Save].
- 6 Tap [Home].
- ⇒ User home screen appears.

Select the method for measuring


- 1 Tap [Measurement].
 - ⇒ The methods list is displayed.
- 2 Tap My first method.
 - ⇒ The work screen of the method **My first method** appears.
 - ⇒ The sample chamber opens automatically.

Placing sample pan

- ► The display now prompts you to load the empty sample pan and tare the balance.
- 1 Place the empty sample pan in the sample pan handler.
- 2 Place the sample pan handler in the sample chamber. Ensure that the tongue of the pan handler fits exactly in the slot of the draft shield element. The sample pan must lie flat in the pan holder.

Note

We advise you to work with the sample pan handler at all times. The pan handler is ergonomic, automatic positioning, safe and provides protection against possible burns due to the hot sample pan.

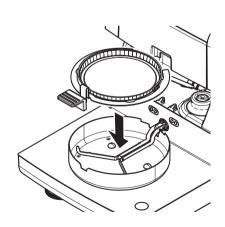


Taring balance

- Tap [->0/T<-].
 - ⇒ The sample chamber closes automatically for taring.
 - ⇒ When taring is completed, the sample chamber opens automatically.

→0/T←

Start measuring

- ► Following taring, the display prompts you to add the sample to the sample pan.
- 1 Place the specimen sample (glass fibre filter) in the sample pan.
- 2 Wet the specimen sample with a few drops of water.
- 3 Tap [Start drying]
 - \Rightarrow The sample camber closes automatically.
 - ⇒ The drying process starts automatically.

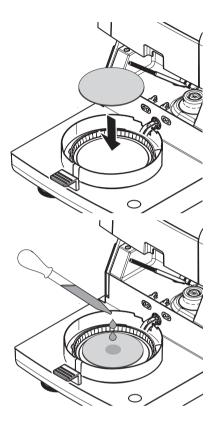

Drying process

You can follow the measurement process on the display, **see** Work Screen (page 21).

- The drying process is continuously displayed graphically.
- The current temperature in the heating module is displayed as well as the elapsed drying time and the current drying value.
- The display shows the selected settings.
- The drying process can be aborted by tapping [Stop drying].

At the end of the drying process you can read the moisture content of your sample on the display. If you have defined the control limits in the method, passed, warned or failed status is shown.

Remove Sample



CAUTION

Danger of burns


Sample, sample pan and sample pan holder may still be hot.

- ▶ The drying process is completed.
- ► The sample chamber is open (opens automatically after the measurement is completed).
- 1 Carefully remove the sample pan handler from the sample chamber.

Note

To remove the sample pan from the handler, lift the pan slightly from below and remove it from the handler.

- 2 To perform a further measurement with the current method, tap [Next Sample].
 - To perform a measurement with a new method, tap [Measurement].
 - To return to the home screen, press [1].

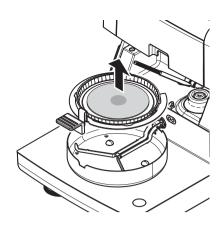

Congratulations!

You have just performed your first measurement with your new Moisture Analyzer. In the following sections you will find more information about the extensive functions of this instrument.

6 Settings

Navigation: Home > Settings

In this section you will find information how to define the general settings of your instrument.

Menu structure

Main menu	Submenu	Further information	
User preferences	Languages	see Languages (page 33)	
	Screen	see Screen (page 33)	
	Password	see Password (page 33)	
System settings	rem settings Regional Settings see Regional Settings		
	Peripherals	see Peripherals (page 34)	
	Default settings	see Default Settings (page 35)	
	Touch screen adjustment	see Touch Screen Adjustment (page 36)	
Application settings	Identifications management	see Identifications Management (page 38)	
	Print and output management	see Print and Output Management (page 38)	
Quality management	Test / Adjustment settings	see Test / Adjustment settings (page 40)	
User management	Groups	see Groups (page 44)	
	Users	see Users (page 46)	
	Account policies	see Account Policies (page 47)	
System and data management	Export / Import	see Export / Import (page 48)	
	Back up / Restore	see Back up / Restore (page 49)	
	Reset	see Reset (page 50)	
	Update	see Update (page 50)	

6.1 User Preferences

Navigation: Home > Settings > User preferences

This menu item defines the basic settings for each user. This settings are stored under the active user profile and apply when working within this profile. When a user profile is called up (login user account), the relevant settings are loaded automatically. **See** User Management (page 44).

You can define the following parameters:

Menu item	Explanation	Further Information
Languages	Defines your preferred display and keyboard lan-	see Languages (page 33)
	guage.	
Screen	Defines the brightness and the color scheme of the	see Screen (page 33)
	display	

6.1.1 Languages

Navigation: Home > Settings > User preferences > Languages

Select your preferred language for operating the terminal.

Factory setting: Generally, the language of the destination country (if available) or English is set.

You can define the following parameters:

Parameter	Explanation	Values
Display lan- guage	Defines user-specific screen language.	English Deutsch Italiano Français Español Português Polski Magyar Český Русский 日本語 中文 한국어/조선말
Keyboard lan- guage	Defines user-specific layout of the Keypad.	English Deutsch Italiano Français Español Português Polski Magyar Český

6.1.2 Screen

Navigation: Home > Settings > User preferences > Screen

You can define the following parameters:

Parameter	Explanation	Values
Brightness	Defines user-specific brightness of the display.	10100 % (80 %)*
Color scheme	Defines user-specific display color scheme.	Blue* Green Orange Grey

^{*} Factory setting

6.1.3 Password

Navigation: Home > Settings > User preferences > Password

In this menu item you can change the user password. It is also possible to create a password for protection or to deactivate the password protection.

You can define the following parameters:

Parameter	Explanation	Values
Password	Change an existing or create a new password. OFF = no pass-	OFF any
	word protection.	

6.2 System Settings

Navigation: Home > Settings > System settings

This menu item defines the basic settings for the system. The system settings apply to the entire system and therefore to all user profiles and applications.

You can define the following parameters:

Menu item	Explanation	Further Information
Regional set- tings	Defines the settings of date & time and keypad country options.	see Regional Settings (page 34)
Peripherals	Defines the configuration of printers and other devices.	see Peripherals (page 34)
Default settings	Defines the default settings of your instrument.	see Default Settings (page 35)
Touch screen adjustment	Adjusts the touch screen.	see Touch Screen Adjustment (page 36)

6.2.1 Regional Settings

Navigation: Home > Settings > System settings > Regional Settings

This menu item defines the settings of date & time.

You can define the following parameters:

Parameter	Explanation	Values
Date	Sets the date of the clock of the instrument.	Day Month Year
Time	Sets the clock time of the instrument.	Hours Minutes
Date format	Defines the date format.	D. MMM YYYY*
	D = Day	MMM D YYYY I
	M = Month	DD.MM.YYYY
	Y = Year	MM/DD/YYYY
		YYYY/MM/DD
		YYYY-MM-DD
Time format	Defines the time format.	24:MM* 12:MM
	M = Minutes	24.MM 12.MM

^{*} Factory setting

6.2.2 Peripherals

Navigation: Home > Settings > System settings > Peripherals

In this menu item you can configure one printer and one host connection.

Remote clients

The MT-SICS host connection (remote client) is enabled by default and only usable with the METTLER TOLEDO specific protocol MT-SICS.

Commands and further information please refer to the Reference Manual MT-SICS Interface Commands for Halogen Moisture Analyzers downloadable from the internet.

<u>www.mt.com/moisture</u>

Printers

No printer is installed by default. To install a printer:

- 1 Tap [**New...**]
 - ⇒ **Peripheral type** list appears.
- 2 Select device.

Note

The only supported printer is RS-P42 Strip Printer.

- 3 Set the parameters.
- 4 To store the settings, tap [Save].
- 5 To activate the interface, tap [ON].
 - ⇒ Status changes from **Disabled** to **Enabled**.

Note

- A printer or host connection can also be deleted.
- For printer settings see Recommended Printer Settings (page 105).
- For defining the printouts see Print and Output Management (page 38).

You can define the following parameters:

Parameter	Explanation	Values
Device Name	Defines a name for the device.	any
Device Port	Defines the device port. Note It is possible to use an USB-to-Serial adapter. If such an adapter is connected, these devices are shown as e.g.SOU1, SOU2 (SOU stands for serial over USB)	COM1*
Baudrate	Defines the speed of the data transmission (date transfer rate / Baud rate).	1200 2400 4800 9600* 19200 38400 57600 115200
DataBits	Defines the number of data bits.	7 8*
Parity	Defines the variant of parity bit.	None* Odd Even
StopBits	Defines the stop bits of the transmitted data.	1 StopBit* 2 StopBits
Flow Control	Defines the type of flow control (Handshake).	None Hardware Xon/Xoff*

^{*} Factory setting

6.2.3 Default Settings

Navigation: Home > Settings > System settings > Default settings

Define the default settings of your instrument.

You can define the following parameters:

Menu item	Explanation	Further Information
Default lan-	Defines a default display and keyboard language.	see Default Languages (page 35)
guages		
Default screen	Defines the default brightness and color scheme.	see Default Screen Settings
settings		(page 36)

6.2.3.1 Default Languages

Navigation: Home > Settings > System settings > Default settings > Default languages

Select the default language for operating the terminal. Factory setting: Generally, the language of the destination country (if available) or English is set.

You can define the following parameters:

Parameter	Explanation	Values
Default display language	Defines default screen language.	English Deutsch Italiano Français Español Português Polski Magyar Český Русский 日本語 中文 한국어/조선말
Default key- board language	Defines default layout of the keyboard.	English Deutsch Italiano Français Español Português Polski Magyar Český

6.2.3.2 Default Screen Settings

Navigation: Home > Settings > System settings > Default settings > Default screen settings

You can define the following parameters:

Parameter	Explanation	Values
Brightness	Defines default brightness of the display.	10100 % (80 %)*
Color scheme	Defines default display color scheme.	Blue* Green Orange Grey

^{*} Factory setting

6.2.4 Touch Screen Adjustment

Navigation: Home > Settings > System settings > Touch screen adjustment

If the alignment of the touch-sensitive areas of the screen correspond not exactly with the position of the buttons e.g. after a software update was performed, then you can solve this error by executing this function.



CAUTION


Do not touch the touch screen with pointed or sharp objects!

This may damage the touch screen.

- ▶ Menu **System settings** appears.
- 1 Choose Touch screen adjustment and tap [Start].
 - ⇒ Adjustment screen appears.
- 2 Touch and hold a stylus on the center of the target until it moves to the next position on the screen.
- 3 Repeat this procedure until a message appears (New calibration settings have been measured).
- 4 Tap the screen to confirm or wait for 30 seconds to cancel (time-out).
 - ⇒ A message box appears.
- 5 To store the setting, tap [OK].

Note

- Tap the center of the target as exact as possible otherwise the process has to be repeated several times.
- Make sure not to touch any other locations on the screen and the fix key areas while doing this adjustment.
- Make sure not touch the screen with your hand.
- Ensure minimum viewing angle.
- The adjustment process cannot be aborted.

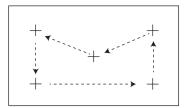

6.3 Applications Settings

Navigation: Home > Settings > Application settings

You can define the following parameters:

Menu item	Explanation	Further Information
Identifications	Defines identifications and their use in order to	see Identifications Management
management	comment measurements.	(page 38).
Print and output	Defines the default settings for the printouts.	see Print and Output Management
management		(page 38).

6.3.1 Identifications Management

Navigation: Home > Settings > Application settings > Identifications management

Identifications (ID) contain descriptive text for individual measurements, which enables samples to be easily allocated to specific tasks or customers. This feature allows you to define identifications in order to comment measurements, such as company id, id badge, id sample. In the ID panel on the work screen, the values (comments) can be entered or changed before measurement, **see** Work Screen (page 21). IDs have an impact on all measurements, regardless of methods. The IDs are included on printouts or can be exported to a PC.

Menu structure

Main menu	Submenu	Further information
Identifications usage	Identifications input	see below
	Clear input after measurement	
Identifications	no submenu	

You can define the following parameters:

Parameter	Explanation	Values
Identifications	Defines the method for entering comments.	OFF* ON
input	ON = activates ID panel on work screen.	Free input Guided input
	Free input = the IDs appear after tapping on the ID panel (work screen) for entries or changes of values (comments).	Outubu IIIpui
	Guided input = each ID is displayed individually in order of definition (ID1 to ID10) before measurement, to make entries or changes of values (comments).	
Clear input after measurement	Defines the ID whose value (comment) is deleted after measurement.	OFF* ON D selection:
Identifications	Defines general identifications in order to comment measurements.	+ -
	+ = New identifications (max. 10 IDs) - = Remove identifications (selectable)	

^{*} Factory setting

6.3.2 Print and Output Management

Navigation: Home > Settings > Application settings > Print and output management

If a printer is connected and activated, results can be automatically printed out. The interim values are printed out in the pre-selected intervals and the final results are logged at the completion of the measurement. The printout contains information to the instrument, time, date as well as the method and results, depending of the protocol. The level of detail depends on the selected type of printout (short, standard) and the activated protocol items (settings or method), **see** Information on Printouts (page 86).

Note

- The definitions in this menu item are the default settings for the printouts. These parameters are displayed at every initialization of a Printout and can be modified again.
- The setting for printout language in this menu item has no impact on test or adjustment printouts, **see** Print and Output Settings (page 42).
- Not all languages are compatible with some printers, and also depending on the settings of the printer, see Recommended Printer Settings (page 105).
- For installing and defining a printer **see** Peripherals (page 34).

Parameter	Explanation	Values
Printout type	Selects a detailed printout or in a short form.	Standard* Short
Printout lan-	Defines the language for general printouts.	English Deutsch
guage	Note No impact on test or adjustment printouts.	Italiano Français Español Português Polski Magyar Český
Print results automatically after measure- ment	Print results automatically after measurement if activated.	(deactivated)* (activated)
Include interme- diate results to printout	Defines the time intervals at which interim results are transferred to the printer.	OFF* 10 s 30 s 1 min 2 min 10 min

^{*} Factory setting

6.4 Quality Management

Navigation: Home > Settings > Quality management

In this menu you can define all settings related to the adjustment and Test of the instrument. Periodic adjustments or tests ensure accurate measurement results.

You can define the following parameters:

Menu item	Explanation	Further Information
Test / Adjust-	Defines the equipment for adjusting and the work-	see Test / Adjustment settings
ment settings	flow for testing or adjusting.	(page 40)

6.4.1 Test / Adjustment settings

Navigation: Home > Settings > Quality management > Test / Adjustment settings

This section describes all the menu options available for defining the parameters relating to the adjustment and testing of your instrument.

For performing test and adjustments see Test/Adjust (page 52).

Menu structure

Main menu	Submenu	Further information
Equipment	Test weight 1	see Equipment (page 40)
	Test weight 2	
	Temperature kit	
Test settings	Weight test - internal	see Test Settings (page 41)
	Weight test - external	
	Temperature test	
Adjustment settings	Weight adjustment - internal	No adjustable settings available
	Weight adjustment - external	see Adjustment Settings (page 42)
	Temperature adjustment	
Print and output settings	Printout language	see Print and Output Settings
	Print results automatically after	(page 42)
	test/adjustment	
FACT	no submenu	see FACT (Fully automatic adjustment) (page 43)

6.4.1.1 Equipment

Navigation: Home > Settings > Quality management > Test / Adjustment settings > Equipment

This menu item defines the equipment needed for weight adjustment, weight test, temperature adjustment or temperature test.

Note

For performing temperature adjustment or temperature test, you need the optional temperature adjustment kit, see Accessories and Spare Parts (page 100).

Test Weight 1 and Test Weight 2

You can define the following parameters:

Parameter	Explanation	Values
Temperature kit	Defines a name that is in an easily recognizable form as an alter-	any
name	native to the ID. A name should be unique and unambiguous.	
Nominal weight	Defines the nominal weight in grams.	30.000200.000 g (200 g)*
Identification	Defines the identification (ID) of the weight. The ID of the weight is provided on the Weight Certificate. The ID may contain your company specific identification number. This ID is stored in the adjustment or test history and shows up on the printouts.	any

^{*} Factory setting

Temperature Adjustment Kit

You can define the following parameters:

Parameter	Explanation	Values
Temperature kit	Defines a name that is in an easily recognizable form as an alter-	any
name	native to the ID. A name should be unique and unambiguous.	
Identification	Defines the identification (ID) of the Temperature kit. The ID is	any
	provided on the Certificate. The ID may contain your company	
	specific identification number. This ID is stored in the adjustment	
	or test history and shows up on the printouts.	

6.4.1.2 Test Settings

Navigation: Home > Settings > Quality management > Test / Adjustment settings > Test settings

Weight test internal

You can define the following parameters:

Parameter	Explanation	Values
Tolerance	Defines the measuring tolerance of the internal weight.	01 %
		(0.001 %)*

Weight test external

You can define the following parameters:

Parameter	Explanation	Values
Test weight	Selects the predefined Test weight . Test weight 1 / Test weight 2 = defined in the menu item Equip-	Test weight 1
	ment	resi weigili 2
Tolerance	Defines the measuring tolerance of the weight.	0.00010.1000 g
		(0.0010 g)*

^{*} Factory setting

Temperature test

Parameter	Explanation	Values
Temperature kit	Shows the name of the Temperature kit which was defined in the	none
	menu item Equipment .	
Temperature 1	Defines the lower test temperature.	OFF ON
	The value is based on the lower adjustment temperature and is	(100 °C)*
	not editable in the temperature test.	
Tolerance	Defines the tolerance of the test temperature.	15 °C
	·	(3 °C)*
Temperature 2	Defines the upper test temperature	OFF ON
	The Value is based on the upper adjustment temperature.	(160 °C)*

Tolerance	Defines the tolerance of the test temperature.	15 °C
		(3 °C)*
Temperature 3	Defines an additional Test temperature.	OFF* ON
(additional)		60200 °C
Tolerance	Defines the tolerance of the test temperature.	15 °C
	·	(3 °C)*

^{*} Factory setting

6.4.1.3 Adjustment Settings

Navigation: Home > Settings > Quality management > Test / Adjustment settings > Adjustment settings

Weight adjustment internal

No editable parameters.

Weight adjustment external

You can define the following parameters:

Parameter	Explanation	Values
Test weight	Selects the predefined Test weight .	Test weight 1
	Test weight 1 / Test weight 2 = defined in the menu item Equip-	Test weight 2
	ment	

Temperature adjustment

You can define the following parameters:

Parameter	Explanation	Values
Temperature kit	Shows the name of the Temperature kit which was defined in the menu item Equipment .	none
Temperature 1	Defines the lower temperature.	80140 °C (100 °C)*
Temperature 2	Defines the upper temperature	140200 °C (160 °C)*

^{*} Factory setting

Note

Most samples are measured in the range of 100 ° C and 160 ° C. This is also the factory setting. The range can be expanded to the range of 80°C and 200°C. The difference between the lower and upper temperature has to be equal or more than 60 °C.

6.4.1.4 Print and Output Settings

Navigation: Home > Settings > Quality management > Test / Adjustment settings > Print and output settings

This menu item allows you to set specific printer functions for test or adjustment printouts.

- The setting for printout language in this menu item has impact on test or adjustment printouts only. For setting the general printout language, see Print and Output Management (page 38).
- Not all languages are compatible with any printer, and also depending on the settings of the printer, see Recommended Printer Settings (page 105).
- For installing and defining a printer **see** Peripherals (page 34).

You can define the following parameters:

Parameter	Explanation	Values
Printout lan- guage	Defines the printout language for test or adjustment printouts.	English Deutsch Italiano Français Español Português Polski Magyar Český
Print results automatically after test/adjust- ment	Activated or deactivated function.	✓ (enabled)* I ☐ (disabled)

^{*} Factory setting

6.4.1.5 FACT (Fully automatic adjustment)

Navigation: Home > Settings > Quality management > Test / Adjustment settings > FACT

This menu item allows you to deactivate or activate fully automatic adjustment function FACT.

Note

- Before performing an external weight adjustment, we recommend you to disable FACT. Otherwise FACT
 overwrites the adjustment.
- This function is not available for approved models that is, **FACT** is always active.

Parameter	Explanation	Values
FACT	Enabled or disabled fully automatic adjustment FACT function.	✓ (enabled)* I☐ (disabled)

^{*} Factory setting

6.5 User Management

Navigation: Home > Settings > User management

The security system of the instrument is based on a "Users and Groups" concept. Every real user of the instrument should be represented by a user account on the instrument which holds particular settings of the respective user. Every user belongs to a particular group and inherits the access rights assigned to that group.

You can define the following parameters:

Menu item	Explanation	Further Information
Groups	Defines groups and their access rights.	see Groups (page 44)
Users	Defines the user accounts.	see Users (page 46)
Account policies	Set default startup user.	see Account Policies (page 47)

6.5.1 Groups

Navigation: Home > Settings > User management > Groups

In this menu you can create, delete, or edit user groups and define their access rights.

Four predefined groups are available. The default group with all access rights is administrator. The administrator group can not be edited or deleted. All other groups can be edited or deleted.

Predefined groups	Default access rights						
	User Pref- erence set- tings	System set- tings	Edit Meth- ods	Run Tests	Run Adjust- ments	Quality manage- ment	User and data management
Administra- tor	V	V	V	V	\checkmark	V	V
Supervisor	V	V	\checkmark	$\overline{\checkmark}$	$\overline{\checkmark}$		
Quality Manager	V	V				V	
Operator	V						

Creating a new group

- ▶ Menu **Groups** selected.
- 1 Tap [**New...**].
- 2 Enter the name of the new group and confirm with [OK]. Name should be unique and unambiguous.
- 3 Edit the properties of the new group.

Deleting a group

- ▶ Menu Groups selected.
- 1 Select the group you wish to delete.
- 2 Tap [Delete]. Confirm the message box.

Note

The Administrator group as well as groups with assigned users cannot be deleted, see Users (page 46).

Editing a group

- ► Menu **Groups** selected.
- 1 Select the group you wish to edit.
- 2 Edit the group parameters. Store the settings with [Save].

You can define the following parameters:

Parameter	Explanation	Values
Group properties	Name and (optional) description of the group. Name should be unique and unambiguous.	any
Access rights	Defines the access rights to menus or operations.	User Preference set- tings System settings Edit Methods Run Tests Run Adjust- ments Quality Man- agement User and data management
Group members	Displays the users assigned to this group (no editing can be done here, the assignment is done when setting up or editing a user).	none
History	Displays the date and time of the last modifications made to this group (no editing here).	none

Influence of access rights

Access rights	Impact on
User Preference settings	User preferences
	Navigation: Home > Settings > User preferences
	See User Preferences (page 33)
System settings	System settings
	Navigation: Home > Settings > System settings
	See System Settings (page 34)
Edit Methods	Method definition
	Navigation: Home > Method definition
	See Method Definition (page 61)
	Application settings
	Navigation: Home > Settings > Application settings
	See Applications Settings (page 38)
Run Tests	Tests
	Navigation: Home > Test/Adjust > Tests
	See Tests (page 57)
Run Adjustments	Adjustments
	Navigation: Home > Test/Adjust > Adjustments
	See Adjustments (page 52)
Quality management	Quality management
	Navigation: Home > Settings > Quality management
	See Quality Management (page 40)
User and data management	User management
	Navigation: Home > Settings > User management
	See User Management (page 44)
	System and data management
	Navigation: Home > Settings > System and data management
	See System- and Data Management (page 48)
	Deleting results
	See Graphical Evaluations of Measurement Results (page 84)

Note

- If password protection is mandatory requirement, the access right User Preference settings must be disabled
- If a user does not have the appropriate access rights, the respective navigation elements are not shown (on Home screen) or are disabled (greyed out).

6.5.2 Users

Navigation: Home > Settings > User management > Users

In this menu you can create, edit, or delete user accounts.

Attention

Important note for the Administrator password.

We recommend that you note down your administrator password and keep it in a safe place. If you forgot the password for the default administrator access, it is not possible to regain access to the moisture analyzer. In this case contact your METTLER TOLEDO representative.

Creating a new user account

Note

The user account name must be unique.

- ▶ Menu **Users** selected.
- 1 Tap [**New...**].
- 2 Enter the name of the new user account and confirm with [OK].
- 3 Edit the properties of the new user account.

Editing an existing user account

- ► Menu **Users** selected.
- 1 Select the user account you wish to edit.
- 2 Select menu User properties.
- 3 Edit the user account parameters.
- 4 To store the settings, tap [Save].

You can define the following parameters:

Parameter	Explanation	Values
User name	Name of the user account. Name should be unique and unambiguous.	any
Full name	Full name of the user (optional).	any
Description	Descriptive text for the user account (optional).	any
Account active	Activate or deactivate the user account. Inactive user accounts are not available for login.	☐ (deactivated)* I ☑ (activated)
Password	Create a new or change an existing password. OFF = no password protection.	OFF* ON any
Group	Assign the user account to a particular group.	available groups

^{*} Factory setting

Note

- To view the date and time of last modification made to a user account, select the respective account and tap [History].
- If password protection is mandatory requirement, the access right User Preference settings must be disabled, see Groups (page 44).

Deleting a user account

- ► Menu **Users** selected.
- 1 Select the user account you wish to delete.
- 2 Tap [**Delete**]. Confirm the message box.

Note

You cannot delete the Administrator account or the account that is currently active.

6.5.3 Account Policies

Navigation: Home > Settings > User management > Account policies

In this menu you can set the default startup user.

You can define the following parameters:

Parameter	Explanation	Values
Default startup user	Select the default user account for login at startup.	OFF ON* available user accounts

^{*} Factory setting

Note

If a non-standard user logs out, the default startup user profile is automatically activated.

6.6 System- and Data Management

Navigation: Home > Settings > System and data management

This menu item allows you to export or import settings and methods to or from an external USB storage device, as well as to restore the system (depending on the user rights).

Menu structure

Main menu	Submenu	Further information
Export / Import	Export settings and methods to	see Export / Import (page 48)
	external storage devices	
	Import settings and methods from	
	external storage device	
Back up / Restore	Back up your system on external	see Back up / Restore (page 49)
	storage device	
	Restore your system from selected	
	back up	
Reset	no submenu	see Reset (page 50)
Update	Update your instrument software	see Update (page 50)

6.6.1 Export / Import

Navigation: Home > Settings > System and data management > Export / Import

This menu item allows you to export or import settings and methods to or from an external storage device. These data can be transferred to other instruments of the same type and can be useful to configure multiple instruments with the same specifications.

Note

- Software Version must be equal or higher.
- Data transfer from HX models to HS models is limited. Features which are not supported by HS models cannot be imported, e.g. Step drying or high resolution.
- Export / import of methods can also be performed from the menu Method Definition.

Export settings and methods to external storage devices...

Parameter	Explanation	Values
File name	Defines the name of the new data file.	any
Location	Finds the location on the storage device.	Browse and select
Export selection	Defines the data for export.	AII* ✓ User manage-
	Note	ment and user prefer- ences
	User management and user preferences contains: User management settings, user preferences	System, application and QM settings
	System, application and QM settings contains: System settings, application settings, quality management settings	✓ Methods

^{*} Factory setting

- ▶ An external storage device is connected e.g. memory stick.
- Settings are made.
- To start, tap [Export].
 To abort, tap [Cancel].

Import settings and methods from external storage device...

You can define the following parameters:

Parameter	Explanation	Values
Import from file	Defines the data to be imported.	Browse and select
Location	Shows the location of the import file.	_
Import selection	Defines the data for import.	AII* ✓ User manage-
		ment and user prefer-
		ences
		and QM settings
		✓ Methods

^{*} Factory setting

- ▶ Import settings and methods from external storage devices is selected.
- ▶ An external storage device is connected e.g. memory stick.
- 1 Tap Browse and select
 - ⇒ **Import from file** appears.
- 2 Select Device and import file.
 - ⇒ Full menu of Import settings and methods from external storage devices appears.
- 3 Tap Import selection and select the data for import if needed.
- 4 To start, tap Import button and follow the instructions in the message window. To abort, tap [Cancel].

6.6.2 Back up / Restore

Navigation: Home > Settings > System and data management > Back up / Restore

This menu item allows you to create a recovery point of the system and to store them on an external storage device. We advise you to create a recovery point of the system regularly. So you are able to restore the system with all data in fault.

Attention

- After restoring, the systemis in backup status. That means, that newer data such as settings, methods or results are lost.
- Restore is only possible on the same instrument as the back up was carried out.

Backup

You can define the following parameters:

Parameter	Explanation	Values
File name	Defines the name of the new data file.	any
Location	Finds the location on the storage device.	Browse and select

Procedure:

- ► An external storage device is connected e.g. memory stick.
- ▶ Back up your system on external storage device... is selected.
- 1 Define the name of the new data file.
- 2 Define the location on the storage device.
- 3 To confirm, tap [OK].
- 4 To start, tap [Backup]. To abort, tap [Cancel].

Restore

You can define the following parameters:

Parameter	Explanation	Values
Restore from file	Finds the location on the storage device and selects the file for	Browse and select
	restoring.	

Procedure:

- ▶ Restore your system from selected back up... is selected.
- ▶ An external storage device is connected e.g. memory stick.
- 1 Tap Restore from file... (Browse and select).
 - ⇒ Storage device appears.
- 2 Select the file for restoring.
- 3 To start, tap [Restore]. To abort, tap [Cancel].
 - ⇒ Confirmation window appears.
- 4 Confirm with [Restore].
- ⇒ The system will restore the data and then perform a restart. Wait until the system has completely started up again.

6.6.3 Reset

Navigation: Home > Settings > System and data management > Reset

This menu item allows you to reset the Instrument to factory settings. Date, time, weight and temperature adjustments/history are not affected.



CAUTION

Data loss after system reset.

After resetting, the instrument is in state as delivered. That means, that all data such as settings, methods or results are lost.

We recommend you, first save the appropriate data on an external storage device, see
 Export / Import (page 48) and Back up / Restore (page 49).

You can define the following parameters:

Parameter	Explanation	Values
Reset	Resets the instrument to factory settings.	none

6.6.4 Update

Navigation: Home > Settings > System and data management > Update

METTLER TOLEDO is continuously improving its instrument firmware (software) for the benefit of customers. So that the customer can benefit quickly and easily from further developments, METTLER TOLEDO makes the latest firmware versions available on the Internet. The firmware made available on the Internet has been developed and tested by Mettler-Toledo AG using processes that meet the guidelines of ISO 9001. Mettler-Toledo AG does not, however, accept liability for consequences that might arise from using the firmware.

Operating Principle

If you register on the following METTLER TOLEDO website, you will find all relevant information and updates for your instrument:

www.mettler-toledo-support.com

Update Procedure



CAUTION

Data loss after system update.

During update the system, the existing data such as settings, methods or results are migrated. It may happen that data are lost or the update fails.

We advise you, first create a recovery point of the system on an external storage device,
 see Back up / Restore (page 49).

For any software update, four steps must be carried out:

Step 1 Backup system data.

See Back up / Restore (page 49), procedure Backup.

Step 2 Execute software update (see below).

Step 3 Restore system data.

See Back up / Restore (page 49), procedure Restore.

Step 4 Adjust the touch screen if necessary (see note).

See Touch Screen Adjustment (page 36).

Execute software update

This menu item allows you to update the Instrument.

You can define the following parameters:

Parameter	Explanation	Values
Update from file	Finds the location on the storage device and selects the file for	Browse and select
	updating.	

- ▶ Back up of the system data is already carried out (Step 1)
- 1 Insert USB stick with software for update.
- 2 Select Home > Settings > System and data management > Update.
 - ⇒ **Update** screen appears.
- 3 Tap **Update from file. Browse and select** the software update package file on the USB stick.
- 4 Wait until the software update is completely finished, i.e. the system returned to the home screen (the instrument might reboot up to twice in between).

Attention

Do not remove the USB stick before the instrument has completely restarted.

5 Restore system data (Step 3).

Note

We recommend to perform a restore. It can be that data are lost.

6 Perform **Touch screen adjustment** if necessary (see note).

Navigation: Home > Settings > System settings > Touch screen adjustment Note

May be necessary if the update also includes the operating system and the alignment of the touch-sensitive areas of the screen correspond not exactly anymore.

7 Test/Adjust

Navigation: Home > Test/Adjust

This function offers you to adjust or test the integrated balance and the heating module. For the relevant settings see Test / Adjustment settings (page 40).

Menu structure

Main menu	Submenu	Further information
Adjustments	Weight adjustment - internal	see Weight Adjustment - Internal
		(page 53)
	Weight adjustment - external	see Weight Adjustment - External
		(page 53)
	Temperature adjustment	see Temperature Adjustment
		(page 54)
	Adjustment Printouts	see Adjustment Printout Examples
		(page 56)
Tests	Weight test - internal	see Weight test - internal (page 57)
	Weight test - external	see Weight test - external (page 57)
	Temperature test	see Temperature test (page 58)
	Test Printouts	see Test Printout Examples
		(page 59)
History	no submenu	see History (page 60)

7.1 Adjustments

Navigation: Home > Test/Adjust > Adjustments

In this menu item you can adjust the balance as well as the heating module of your instrument. For relevant settings **see** Test / Adjustment settings (page 40).

Attention

- To obtain accurate results, the balance must be adjusted at the point of use under measuring condition in order to match the gravitational acceleration at its location. The instrument must be connected to the power supply for approximately 60 minutes in order to reach operating temperature before adjusting. Adjusting is necessary:
 - · before the instrument is used for the first time.
 - after a change of location.
- We advice you to wait at least 30 minutes after a drying operation (or previous heating module adjustment) before performing the adjustment.

Fully automatic adjustment FACT

This instrument is equipped with fully automatic adjustment with two internal weights called Fully Automatic Calibration Technology (**FACT**).

The balance adjusts itself automatically:

- after the warm-up phase after connection to the power supply.
- when a change in the ambient conditions (temperature) could lead to a noticeable deviation in the measurement.
- on a fixed time interval

FACT is activated by default and can be switched off as required e.g. when performing an external adjustment, see FACT (Fully automatic adjustment) (page 43).

Navigation: Home > Settings > Quality management > Test / Adjustment settings > FACT

Note

With approved models **FACT** cannot be switched off.

7.1.1 Weight Adjustment - Internal

Navigation: Home > Test/Adjust > Adjustments > Weight adjustment - internal

This function allows you to adjust the balance with the built-in weights. The internal adjustment is equivalent to the FACT adjustment.

Perform this function as follows:

- 1 Tap the menu item Weight adjustment internal.
 - ⇒ The adjustment screen appears.
- 2 Clear pan and tap [Start].
 - ⇒ The balance adjusts itself automatically.
- ⇒ The instrument indicates when the adjustment process is done. The balance is adjusted and ready to use for further measurements. The display shows the adjustment results. The results can be printed out by tapping [□].

7.1.2 Weight Adjustment - External

Navigation: Home > Test/Adjust > Adjustments > Weight adjustment - external

This function allows you to adjust the balance using an external weight. We recommend to use certified weights so that full traceability is guaranteed.

Note

We recommend you to disable FACT. Otherwise FACT overwrites the adjustment.

Perform this function as follows:

- ▶ FACT is off.
- ► Test weight and Adjustment settings are defined. see Test / Adjustment settings (page 40)
- 1 Have required test weight ready.
- 2 Tap the item Weight adjustment external.
 - ⇒ The adjustment screen appears.
- 3 Clear sample pan holder and tap [Start].

Note

We recommend not to use a pan.

- ⇒ The required value flashes and prompts you to load the weight, e.g. "100.00 g".
- 4 Load the required test weight in the center of sample pan holder.
 - ⇒ The balance indicates the adjustment process, "———" appears.
- 5 When the display flashes "000.00" remove the test weight.
- ⇒ The balance is adjusted and ready to use for further measurements. The display shows the adjustment results. The results can be printed out by tapping [♣]

Note

- After a timeout the error message appears Weight out of range. The adjustment has been aborted The adjustment was not executed.
- 1 Confirm with [**OK**].
- 2 Remove the test weight for unloading.
- 3 Repeat the adjustment.

7.1.3 Temperature Adjustment

Navigation: Home > Test/Adjust > Adjustments > Temperature adjustment

This function allows you to adjust the temperature control of the heating module. You need the optional **Temperature kit** for performing this function, **see** Accessories and Spare Parts (page 100). To learn when an adjustment of the heating module is necessary, **see** Notes on Adjustment of Balance and Heating Module (page 104). We advise you to wait at least 30 minutes after a drying operation (or previous heating module adjustment) before performing the adjustment.

Note

- The duration of the temperature adjustment is 30 minutes. (15 minutes for each measuring temperature).
- The sample chamber must be less than 50 °C before a further adjustment or test can be done.
- It is also possible to perform a test first, and then to transform the test into an adjustment if required. Only possible for tests conducted with 2 temperatures, **see** Temperature test (page 58).
- For temperature adjustment settings see Adjustment Settings (page 42).



CAUTION


Danger of burns

Sample chamber may still be hot.


- Cool down instrument before setting up.
- ▶ Adjustment settings are defined, **see** Adjustment Settings (page 42).
- Tap menu item Temperature adjustment.
 - ⇒ Work screen **Temperature adjustment** appears.

Setting up Drying Unit

- ► Have required **Temperature kit** ready.
- Instrument is on
- ▶ Sample chamber is open.
- ▶ Instrument is cooled down.
- 1 Remove sample pan handler (with pan) (1).
- 2 Remove sample pan holder (2).
- 3 Contact areas (3) should be clean. Clean them if needed.

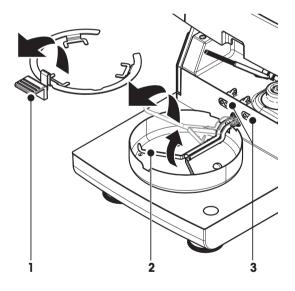

Insert Temperature kit (4).

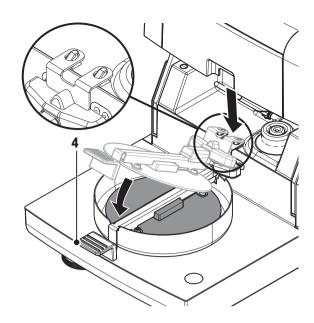

Start adjustment

- ► Temperature kit is inserted.
- 1 Tap Start.

Note

The adjustment process takes 30 minutes.

- ⇒ The instrument starts the adjustment process. The screen shows the current temperature and the remaining time.
- ⇒ The instrument heats up to the target temperature 1 for determining the predefined lower temperature.
- ⇒ After 15 minutes, the instrument adjusts the lower temperature and continues with the process.
- ⇒ The instrument heats up to the target temperature 2 for determining the predefined upper temperature.
- ⇒ After 15 minutes, the instrument adjusts the upper temperature and concludes the adjustment.
- 2 To print out the results, tap [=].
 - To display or print out the results of previous adjustments, tap [History].
 - To return to the home screen, tap [1].
- 3 After cooling down, remove **Temperature kit** and set up drying unit for measurement.

See History (page 60) and Setting up Drying Unit (page 25)

Note

Your calibrated temperature adjustment kit can be re-calibrated. Please contact your local METTLER TOLEDO representative for more information.

7.1.4 Adjustment Printout Examples

Weight adjustment internal

Weight adjustment external

-WEIGHT ADJUST EXTERNAL-
14.10.2011 11:51
METTLER TOLEDO Halogen MoistureAnalyzer
Type HX204 SNR(Drying Unit) 2345 SNR(Terminal) 6788 SW(Drying Un 0.60.09 SW(Terminal) 4.20628
User name Administrator
Weight ID ECW-100/1
Nominal Weight 100.000 g
Cell temperature 25.20 °C
Drying Unit leveled Adjustment done
Signature
END

Temperature adjustment

-TEMPERATURE ADJUSTMENT-
14.10.2011 11:51
METTLER TOLEDO Halogen MoistureAnalyzer
Type HX204 SNR(Drying Unit) 2345 SNR(Terminal) 6788 SW(Drying Unit) 0.60.09 SW(Terminal) 4.20628
User name Administrator
Temp Kit ID MT-414/A Temp1 target 100 °C Temp1 actual 97 °C Temp2 target 160 °C Temp2 actual 162 °C Adjustment done
Signature
END

7.2 Tests

Navigation: Home > Test/Adjust > Tests

In this menu item you can check the adjustment of the balance as well the temperature of the heating module. For relevant settings **see** Test / Adjustment settings (page 40).

7.2.1 Weight test - internal

Navigation: Home > Test/Adjust > Tests > Weight test - internal

This function allows you to test the balance with the built-in weight. Perform this function as follows:

- ➤ Test settings are defined (Factory setting of the tolerance is 0.0010 %) see Test / Adjustment settings (page 40)
- 1 Open sample chamber
- 2 Tap the item Weight test internal.
 - ⇒ The test screen appears.
- 3 Clear pan and tap [Start].
 - \Rightarrow The balance is tested automatically.
- ⇒ When the test process is done, the display shows the test results and indicates whether the test is passed or failed according to defined tolerances.

The test results can be printed out by tapping [].

7.2.2 Weight test - external

Navigation: Home > Test/Adjust > Tests > Weight test - external

This function allows you to test the balance using an external weight. Perform this function as follows:

- Test weight and test settings are defined.
 see Test / Adjustment settings (page 40)
- 1 Have required test weight ready.
- 2 Open sample chamber.
- 3 Tap the item Weight test external.
 - \Rightarrow The test screen appears.
- 4 Clear sample pan holder and tap [Start].

Note

We recommend not to use a pan.

- ⇒ The required value flashes and prompts you to load the weight.
- 5 Load the required test weight.
 - ⇒ The balance indicates the test process, "———" appears.
- 6 When the display flashes "000.00" remove the test weight.
- ⇒ When the test process is done, the display shows the test results and indicates whether the test is passed or failed according to defined tolerances. The results can be printed out by tapping [◄].

Timeout

- ▶ After a timeout of 75 seconds the error message appears **Weight out of range. The test has been aborted.**The test was not executed.
- 1 Confirm with [**OK**].
- 2 Repeat the test.

7.2.3 Temperature test

Navigation: Home > Test/Adjust > Tests > Temperature test

This function allows you to test the temperature control of the heating module. You need the optional Temperature Kit for performing this function, **see** Accessories and Spare Parts (page 100). To learn when an test of the heating module is necessary, **see** Notes on Adjustment of Balance and Heating Module (page 104). We advise you to wait at least 30 minutes after a drying operation (or previous heating module adjustment) before performing a test.

Note

- The duration of the temperature test is 15 minutes for each measuring temperature.
- The sample chamber must be cooled down below 50 °C before a further adjustment or test can be done.
- It is possible to perform a test and then subsequently to transform the test into an adjustment if required. This is only possible for tests conducted with temperatures 1 and 2.
- Temperature test settings see Test Settings (page 41).

Setting up the Drying Unit for temperature test

Same procedure as for temperature adjustment, see Temperature Adjustment (page 54).



CAUTION

Danger of burns

Sample chamber may still be hot.

Cool down instrument before setting up.

Performing temperature test

- ► Test settings are defined, see Test Settings (page 41).
- ▶ Setting up the drying unit for temperature testing has been carried out.
- 1 Tap menu item Temperature test.
 - ⇒ Work screen **Temperature test** appears.
- 2 Tap [Start].

Note

The testing process takes 15 minutes for each temperature (depending on the settings).

- ⇒ The instrument starts the testing process. The screen shows the current temperature and remaining time.
- ⇒ The instrument is heating up to the target temperature 1 for determining the predefined lower temperature.
- After 15 minutes, the instrument displays and stores the measured lower temperature and continues with the process.
- ⇒ The instrument is heating up to the target temperature 2 for determining the predefined upper temperature.
- ⇒ After 15 minutes, the instrument displays and stores the measured upper temperature.
- ⇒ If temperature 3 is defined, the procedure has an additional step.
- ⇒ The test is done and indicates whether the test has passed or failed according to defined tolerances.
- 3 Decide if an adjustment is needed.
- 4 To return to the menu **Tests**, tap [**Tests**].
 - To print out the results, tap [].
 - To display or print the results at any time, tap [History].
 - To return to the home screen, press [1].
- 5 After cooling down, remove temperature adjustment kit and set up drying unit for measurement, **see** Setting up Drying Unit (page 25).

7.2.4 Test Printout Examples

Weight test internal

Weight test external

WEIGHT TEST EXTERNAL
14.10.2011 11:51
METTLER TOLEDO Halogen MoistureAnalyzer
Type HX204 SNR(Drying Unit) 2345 SNR(Terminal) 6788 SW(Drying Unit) 0.60.09 SW(Terminal) 4.20628
User name Administrator
Weight ID ECW-100/1
Nominal Weight
100.000 g Actual 100.001 g Difference 0.001 g Tolerance +/- 0.002 g
Cell temperature
23.83 °C Drying Unit leveled Test passed
Signature
END

Temperature test

TEMPERATURE TEST
14.10.2011 11:51
METTLER TOLEDO Halogen MoistureAnalyzer
narogen norbeatermaryzer
Type HX204
SNR (Drying Unit) 2345
SNR (Terminal) 6788
SW(Drying Unit) 0.60.09
SW(Terminal) 4.20628
User name Administrator
Temp Kit ID MT-414/A
Temp1 target 100 °C
Temp1 actual 99 °C
Temp1 tolerance
+/- 2 °C
Temp2 target 160 °C
Temp2 actual 159 °C
Temp2 tolerance
+/- 2 °C
Test passed
Signature
END

7.3 History

Navigation: Home > Test/Adjust > History

The instrument always records all test or adjustment operations that have been made and saves them in a memory which is specially protected against power failure (limited to the last 50 records per history selection).

Note

Aborted adjustments or tests will not be saved.

When the History is selected, a list appears with the history selection as follows:

- Weight test history
- Weight adjustment history
- Temperature test history
- Temperature adjustment history

The following information is displayed:

Weight test history

Weight test external		Weight test internal	
Calibration weight identification		Nominal weight	100 %
Nominal weight	g	Actual weight	%
Actual weight	g	Difference	%
Difference	g	Tolerance	%
Tolerance	g	Temperature	°C
Temperature	°C	Leveling information	
Leveling information	·		

Weight adjustments history

Weight adjustment external		Weight adjustment internal	
Calibration weight name		Temperature	°C
Calibration weight identification		Leveling information	
Nominal weight	g		
Temperature	°C		
Leveling information			

Temperature adjustment/test history

Temperature adjustment		Temperature test	
Adjustment kit identification		Adjustment kit identification	
Temperature 1	target - actual	Temperature 1	target - actual
Temperature 2	target - actual	Temperature 2	target - actual
		Free temperature	target - actual
		Tolerances for each temperature	°C

The selected individual results can be printed out by tapping [].

8 Method Definition

Navigation: Home > Method definition

What is a method?

Methods simplify and accelerate your daily work. A method contains all settings for measuring the moisture content of a particular sample (substance). It can be recalled by tapping a measurement button or a shortcut and your instrument immediately operates with the corresponding settings.

The **optimal setting of parameters and the drying time** depends on the type and size of the sample and the desired accuracy of the measurement result. The exact parameters can only be determined experimentally, **see** How to Obtain Best Results (page 104). Your Instrument helps you to determine the settings.

Method definition

This function offers you to define new methods and to change, delete, export or import existing methods depending on the user rights. It allows you to store up to 300 individual methods.

All parameters using in a drying method, can be defined under this menu item.

Note

- In most cases it is sufficient to set Main measurement parameters.
- You will find more information about defining methods in the delivered application brochure «Guide to Moisture Analysis» or see How to Obtain Best Results (page 104).

Prerequisite

The parameter menu appears only if a method already exists and is selected in the method definition dialog or create a new method.

Menu structure

Main menu	Submenu	Further information
Main measurement parame-	Drying program	see Settings of Drying Program
ters	(including temperature and switch-off criteria)	(page 63)
	Display mode	see Settings of Display Mode (page 69)
	Start weight	see Settings of Start Weight (page 71)
Result and value handling	Control limits	see Control limits (page 72)
	Resolution	see Resolution (page 73)
Workflow handling	Starting mode	see Starting Mode (page 75)
	Preheating	see Preheating (page 75).
General method properties	Method name	see Method name (page 77)

Defining a new Method

Note

A method name can only exist once.

- Method definition is selected.
- 1 Tap [New...].
 - \Rightarrow The keypad appears.
- 2 Enter a name for the new method. Name should be unique and unambiguous.1 to 30 characters are possible (including spaces).
- 3 Confirm with [OK].
 - ⇒ The parameter menu for the new method appears.
- 4 Set the desired parameters e.g. Main measurement parameters.
- 5 To store the method, tap [Save].

Editing an existing method

- ▶ Method definition is selected.
- 1 Tap the method in the list you want to edit.
 - ⇒ The parameter menu for the selected method appears.
- 2 Edit the desired parameters e.g. Main measurement parameters.
- 3 To store the method, tap [Save].

Copying an existing method

- Method definition is selected.
- 1 Tap the method in the list you want to copy.
 - ⇒ The parameter menu of the desired method appears.
- 2 Tap [Save as...].
 - \Rightarrow The keypad appears.
- 3 Enter a new name for the copied method. Name should be unique and unambiguous.1 to 30 characters are possible.
- 4 To store the method, tap [Save].

Removing an existing method

Note

All results of this method are also removed.

- ▶ Method definition is selected.
- 1 Tap the method in the list you want to remove.
 - ⇒ The parameter menu of the desired method appears.
- 2 Tap [Delete].
 - ⇒ A message box appears.
- 3 Confirm with [Delete].
- ⇒ The desired method is removed.

Note

A possible shortcut to this method is not removed.

For removing shortcuts see Using Shortcuts (page 82)

8.1 Main Measuring Parameters

Navigation: Home > Method Definition > Method name > Main measurement parameters

These parameters are relevant for measuring and must be determined for all samples. Most samples can be determined with these parameters.

You can define the following parameters:

Menu item	Explanation	Further Information
Drying program	Defines the drying program which is best suited for the specific sample. (including temperature and switch-off criteria)	see Setting the Drying Program (page 63)
Display mode	Defines the type of value for displaying and printing.	see Settings of Display Mode (page 69)
Start weight	Defines a start weight so that the samples always have approximately the same weight.	see Settings of Start Weight (page 71)

8.1.1 Settings of Drying Program

Navigation: Home > Method Definition > Method name > Main measurement parameters > Drying program

This function offers you different predefined drying programs for optimum matching of the drying characteristics to the sample used. Furthermore, the drying temperature, switch-off criterion as well as the start weight can be set. Carry out a test measurement for determining the appropriate switch-off criterion, if you do not know the behavior of the sample. For carry out a test measurement **see** Test a Method (page 78).

Note

You will find more information about defining methods in the delivered application brochure «Guide to Moisture Analysis».

You can define the following parameters:

Parameter	Explanation	Values
Drying program	Defines the drying program which is best suited for the specific	Standard* Rapid
	sample.	Gentle Steps

^{*} Factory setting

Further information:

- Standard drying see Settings of Standard drying (page 63)
- Rapid drying see Settings of Rapid drying (page 64)
- **Gentle** drying **see** Settings of Gentle drying (page 65)
- **Steps** drying **see** Settings of Step drying (page 66)

8.1.1.1 Settings of Standard drying



Standard drying

This drying program is set in the factory and is suitable for most samples. The sample is heated to the drying temperature.

Factory setting

Drying temperature = $105 \, ^{\circ}\text{C}$, Switch-off criterion $3 = 1 \, \text{mg} / 50 \, \text{s}$

You can define the following parameters:

Parameter	Explanation	Values
Drying tempera- ture	Defines the drying temperature.	40230 °C
Switch-off crite- rion	Defines the criterion when the instrument should end the drying.	1 (1 mg / 10 s) 2 (1 mg / 20 s) 3 (1 mg / 50 s)* 4 (1 mg / 90 s) 5 (1 mg / 140 s) Timed Free (mg / s) Free (% / s)

Setting the Drying Temperature

Navigation: Method Definition > Name of method > Main measurement parameters > Drying temperature In this menu item you can set the final drying temperature.

Note:

The admissible input range for the drying temperature is indicated.



- 1 Start of drying
- 2 Final temperature

Setting Switch-off criterion

see Settings of Switch-off Criterion (page 67)

8.1.1.2 Settings of Rapid drying



Rapid drying

This drying program is primarily suitable **for samples with a moisture content over 30%**. Following the start, the selected temperature is exceeded by 40% for 3 minutes (however, is possible up to maximum 230°C) to compensate for the cooling due to vaporization and accelerate the drying process. The drying temperature is then lowered to the set value and maintained.

Factory setting

Drying temperature = $105 \, ^{\circ}\text{C}$, **Switch-off criterion** $3 = 1 \, \text{mg} / 50 \, \text{s}$

Parameter	Explanation	Values		
Drying tempera-	Drying tempera- Defines the drying temperature.			
ture				
Switch-off crite- rion	Defines the criterion when the instrument should end the drying.	1 (1 mg / 10 s) 2 (1 mg / 20 s) 3 (1 mg / 50 s)* 4 (1 mg / 90 s) 5 (1 mg / 140 s) Timed Free (mg / s) Free (% / s)		

^{*} Factory setting


Setting the Drying Temperature

Navigation: Method Definition > Name of method > Main measurement parameters > Drying temperature

In this menu item you can set the final drying temperature.

Note:

The admissible input range for the drying temperature is indicated.

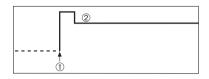

- 1 Start of drying
- 2 Final temperature

Setting Switch-off criterion

see Settings of Switch-off Criterion (page 67)

8.1.1.3 Settings of Gentle drying



Gentle drying

This drying program is suitable for the gentle drying of **substances which tend to form a skin** (e.g. substances containing sugar or volatile substances). With this program the temperature is continuously increased and attains the selected drying temperature **only after the elapse** of the so-called ramp time. The «ramp», i.e. the time that should elapse between the start of drying and attainment of the final temperature can be pre-set.

Factory setting

Drying temperature = 105 °C, Ramp time = 3:00 min, Switch-off criterion 3 = 1 mg / 50 s

You can define the following parameters:


Parameter	Explanation	Values
Drying tempera-	Defines the drying temperature.	40230 °C
ture		
Ramp time	Defines the time in minutes that should elapse between the start of drying and attainment of the final temperature for Gentle drying.	0480 min
Switch-off crite- rion	Defines the criterion when the instrument should end the drying.	1 (1 mg / 10 s) 2 (1 mg / 20 s) 3 (1 mg / 50 s)* 4 (1 mg / 90 s) 5 (1 mg / 140 s) Timed Free (mg / s) Free (% / s)

^{*} Factory setting

Setting the Drying Temperature

Navigation: Method Definition > Method name > Main measurement parameters > Drying programGentleDrying temperature

In this menu item you can set the desire final drying temperature.

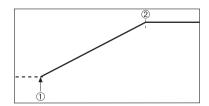


- 1 Start of drying
- 2 Final temperature

Setting the ramp time

Navigation: Method Definition > Method name > Main measurement parameters > Drying programGentle > Ramp time

In this menu item you can define the **Ramp time**, i.e. the time that should elapse between the start of drying and attainment of the final temperature. The ramp is started once the temperature reaches 50 °C.

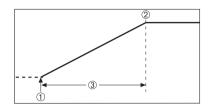

- 1 Start of drying
- 2 Final temperature
- 3 Ramp time

Setting Switch-off criterion

see Settings of Switch-off Criterion (page 67)

8.1.1.4 Settings of Step drying



Step drying

This drying program is suitable for the drying of substances composed of several components which vaporize at different temperatures (e.g. ethereal oils). With this program drying is performed stepwise, i.e. the sample is preheated to a particular temperature (1st step) and kept at this temperature for the selected switch-off criterion e.g. time or weight loss per unit of time. The temperature can be increased for the next step, lowered or remain the same. At the last step, the switch-off criterion terminates the measurement. Maximum 5 steps are selectable.

Step drying can also be used for faster determining of samples with high moisture content. In this case the first step should be set with a higher temperature than for the last step. This reduces evaporation.

Tip for the measurement of plastic granules:

Setting of switch-off criterion "timed" for the first step (possibly same temperature setting as for step 2). The larger sample (e.g. 30 g) has to be heated first before it gives off the small amount of moisture. This prevents that the measurement is terminated prematurely. Setting of switch-off criterion "5" for the second step which terminates the measurement automatically.

Factory setting

Step 1, **Temperature** = $50 \, ^{\circ}$ C, **Switch-off criterion** = $5:00 \, \text{min}$


Step 2, **Temperature** = $105 \, ^{\circ}$ C, **Switch-off criterion** 3 = $1 \, \text{mg} / 50 \, \text{s}$

Parameter	Explanation	Values
Drying steps	ing steps Defines the number of steps for Step drying. This dialog includes the setting of temperature and switch-off criterion for each step.	
Temperature	Defines the drying temperature for each Step.	50230 °C
Switch-off crite- rion	Defines the criterion when the instrument should end the drying or should start the next step.	1 (1 mg / 10 s) 2 (1 mg / 20 s) 3 (1 mg / 50 s)* 4 (1 mg / 90 s) 5 (1 mg / 140 s) Timed Free (mg / s) Free (% / s)

^{*} Factory setting

Drying steps

Navigation: Method Definition > Method name > Main measurement parameters > Drying programSteps > Drying steps

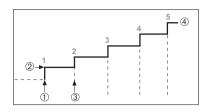

- 1 Start of drying
- 2 Step temperature
- 3 Next step, Switch-off criterion
- 4 Last step

Setting Switch-off criterion

see Settings of Switch-off Criterion (page 67)

8.1.1.5 Settings of Switch-off Criterion

Navigation: Home > Method Definition > Method name > Main measurement parameters > Drying program

This function offers you different switch-off criteria. A switch-off criterion defines when the instrument should end the drying or should start the next step ("Step drying"). Switch-off criteria ensure that measurements are finished in the same condition (weight loss per time) every time, ensuring repeatable measurements.

Note

Accurate results require same start weights, **see** Settings of Start Weight (page 71).

You can define the following parameters:

Parameter	Explanation	Values
Switch-off crite- rion	Defines the criterion when the instrument should end the drying.	1 (1 mg / 10 s) 2 (1 mg / 20 s) 3 (1 mg / 50 s)* 4 (1 mg / 90 s) 5 (1 mg / 140 s) Timed Free (mg / s) Free (% / s)

^{*} Factory setting

- Weight loss per unit of time (5 preprogrammed settings)
- Free switch-off criterion (2 different settings)
- Timed switch-off

Weight loss per unit of time

This switch-off is based on a weight loss per unit of time. As soon as the mean weight loss is less that a preset value during a specified time, the instrument considers drying as complete and automatically discontinues the measurement process.

Note


The switch-off criterion is inactive during the first 30 sec.

The following 5 preprogrammed settings are available

Switch-off criterion	∆g in mg	Δ t in seconds	Description
1	1 mg	10 s	This setting is suitable for fast measurements to determine a trend.
2	1 mg	20 s	This setting is suitable for quick drying samples.
3	1 mg	50 s	This is the factory setting . It is suitable for most kind of samples.

Switch-off criterion	Δg in mg	Δ t in seconds	Description
4	1 mg	90 s	This setting is suitable for samples which dry moderately quickly or for samples with higher accuracy requirements.
5	1 mg	140 s	This setting is suitable for samples which dry very slowly (trapped moisture, skin formation) or for samples with very low moisture content e.g. plastics. Not suitable for very temperature sensitive samples.

The graph below exemplifies the mode of operation of the switch-off operation (not to scale).

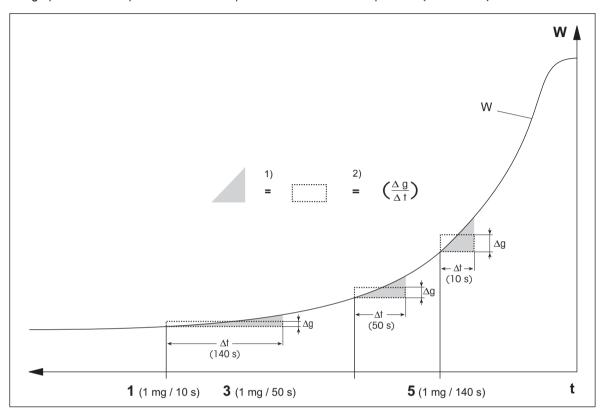

t = time

W = weight of sample

- 1, 3, 5 = switch-off criteria shown as an example
- 1) = equal area
- 2) = mean weight loss per unit of time

Free switch-off criterion

The free switch-off criterion is based on a user-defined mean weight loss per unit of time.

The following 2 settings are available:

- **Switch-off criterion** > **Free (mg / s)...** (weight loss per unit of time)
- **Switch-off criterion** > **Free** (% / s)... (weight loss in percent per unit of time)

Timed switch-off

With this switch-off criterion the measurement lasts until the present drying time has elapsed. The display provides you with continuous information in the drying time.

Performing Test measurements

For performing test measurements to determine a correct switch-off criterion, see Test a Method (page 78).

8.1.2 Settings of Display Mode

Navigation: Home > Method Definition > Method name > Main measurement parameters > Display mode

With this function you can select the desired type of result display. You also define what type of values are printed on the records.

You can define the following parameters:

Parameter	Explanation	Values
Display mode	Defines the type of value for displaying and printing.	%MC* %DC %AM %AD g g/kg MC g/kg DC


- * Factory setting
- %MC Moisture Content
- %DC Dry Content
- %AM ATRO Moisture Content
- %AD ATRO Dry Content (Wet weight)
- **g** Weight in Grams
- g/kg MC Moisture Content
- g/kg DC Dry Content

Detailed information:

%MC - Moisture Content

The moisture content of the sample is displayed (and printed out) as a percentage of the wet weight (WW = initial weight = 100 %). This is the **factory setting**.

During the measuremnet the value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "%MC" (Moisture Content, e.g. 11.35 %MC) also for the printed results.

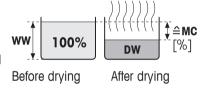

$$MC = \frac{WW - DW}{WW} \cdot 100 \%$$

MC = Moisture Content [0...100 %]


WW = wet weight

DW = dry weight

%DC - Dry Content

The dry content of the sample is displayed (and printed out) as a percentage of the wet weight (WW = initial weight = 100 %).

During the measuremnet the value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "%DC" (Dry Content, e.g. 88.65 %DC) also for the printed results.

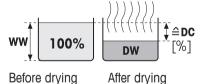

$$DC = \frac{DW}{WW} \cdot 100 \%$$


DC = dry content [100...0 %]

WW = wet weight

DW = dry weight

%AM - ATRO Moisture Content 1)

The moisture content of the sample is displayed (and printed out) as a percentage of the dry weight (DW = final weight = 100 %)

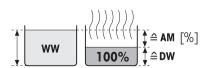

During the measuremnet the value is constantly displayed in percent and Before drying as a graphical drying curve. The measured value is marked by "%AM" (ATRO Moisture Content, e.g. 255.33 %AM) also for the printed results.

$$AM = \frac{WW - DW}{DW} \cdot 100 \%$$

AM = ATRO moisture content [0...1000 %]


WW = wet weight

DW = dry weight

%AD - ATRO Dry Content (Wet weight)1)

The wet weight of the sample is displayed (and printed out) as a percentage of the dry weight (DW = final weight = 100 %)

During the measuremnet the value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "%AD" (ATRO Dry Content, e.g. 312.56 %AD) also for the printed results..

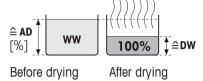

$$AD = \frac{WW}{DW} \cdot 100 \%$$

AD = ATRO dry content [100...1000 %]

WW = wet weight

DW = dry weight

1) Comment on the ATRO display mode

If the current measured value in the ATRO display mode is greater or less than the predefined limit value (i.e. greater than 999.99 %AD or less than –999.99 %AM), the ATRO result values are limited to 999.99%.

g - Weight in Grams


The weight of the sample is displayed (and printed out) in grams. With this setting, the Moisture Analyzer is used as a precision balance.

During the measuremnet current weight is constantly displayed in grams.

g/kg MC - Moisture Content

The moisture content of the sample is displayed (and printed out)in g/kg of the wet weight (WW = initial weight = 1000 g/kg).

During the measuremnet the current measured value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "g/kg MC" (Moisture Content, e.g. 11.35 g/kg MC) also for the printed results.

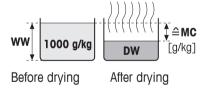

$$MC = \frac{WW - DW}{WW} \cdot 1000 \text{ g/kg}$$


MC = Moisture Content [0...1000 g/kg]

WW = wet weight

DW = dry weight

g/kg DC - Dry Content

The dry content of the sample is displayed (and printed out) in g/kg of the wet weight (WW = initial weight = 1000 g/kg).

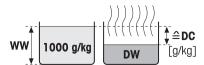

During the measuremnet the current measured value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "g/kg DC" (Dry Content, e.g. 88.65 g/kg DC) also for the printed results.

Before drying After drying

$$DC = \frac{DW}{WW} \cdot 1000 \text{ g/kg}$$

DC = dry content [1000...0 g/kg] WW = wet weight DW = dry weight

8.1.3 Settings of Start Weight

Navigation: Home > Method Definition > Method name > Main measurement parameters > Start weight

The start weight affects both the duration of the measurement as well as the accuracy of the results. A small weight results in a shorter duration of the measurement but will decrease the accuracy of the result. With this function you can define a start weight so that the samples are approximately of equal weight thus improving the repeatability of the measurements. The weighing-in aid assists you in weighing-in the sample. At most of the samples the target weight is in the range of 2–5 g (plastics 30 g). We recommend to cover the whole surface of the sample pan with a thin uniform layer of sample.

Parameter	Explanation	Values
Start weight	Defines a start weight so that the samples always have approxi-	OFF* ON
	mately the same weight .	(0.100200.000 g)
Start weight tol-	Defines the tolerance of the Start weight .	125 %
erance	10 %, recommended for sample weights up to 5 g.	(10 %)*
Weigh-in aid	Defines monitoring of target weight and tolerance.	Passive* Active
	Passive = tolerance is displayed. Active = tolerance is monitored. If the startweight is out of tolerance, the measurement cannot be started.	

^{*} Factory setting

8.2 Result and Values Handling

Navigation: Home > Method Definition > Method name > Result and value handling

You can define the following parameters:

Menu item	Explanation	Further Information
Control limits	Defines the acceptable range for measuring results in the unit of the selected display mode.	see Control limits (page 72)
Resolution	Defines the readability of the built-in analytical balance.	see Resolution (page 73)

8.2.1 Control limits


Navigation: Home > Method Definition > Method name > Result and value handling > Control limits

With this function you can set warning and intervention limits for the quality and process monitoring. You can also set limits for distinguish passed or failed.

Parameter	Explanation	Values
Control limits	Defines the acceptable range for the measuring results in the unit of the selected display mode.	OFF* ON
	OFF = no limits applied.	
T1+ (Upper warning limit)	Defines the upper warning limit in the selected unit of the display mode. Measurement passed with warning (!). Value range in ATRO mode (%AD, %AM) up to 1000.	OFF* 0.01100.0 (1000)
T1- (Lower warning limit)	Defines the lower warning limit in the unit of the selected display mode. Measurement passed with warning (!). Value range in ATRO mode (%AD, %AM) up to 1000.	OFF* 0.01100.0 (1000)
T2+ (Upper intervention limit)	Defines the upper intervention limit in the unit of the selected display mode. Measurement failed (X). Value range in ATRO mode (%AD, %AM) up to 1000.	OFF* 0.01100.0 (1000)
T2- (Lower intervention limit)	Defines the lower intervention limit in the unit of the selected display mode. Measurement failed (X). Value range in ATRO mode (%AD, %AM) up to 1000.	OFF* 0.01100.0 (1000)

^{*} Factory setting

Setting the limits

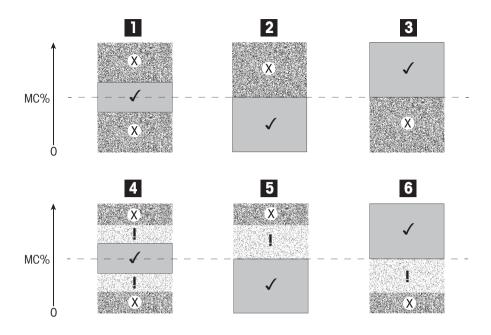

Limits		1	2	3	4	5	6
T2+ (Upper intervention limit)	(X)			Off			Off
T1+ (Upper warning limit)	(!)	Off	Off	Off	0	0	Off
T1- (Lower warning limit)	(!)	Off	Off	Off	0	Off	0
T2- (Lower intervention limit)	(X)		Off			Off	

- (1) Measurement is within the warning limits: passed (displayed in green)
- (!) Measurement is between the warning limits and the intervention limits: passed with warning (displayed in yellow)
- (X) Measurement is outside the intervention limit: failed (displayed in red)

Control limits is off = no limits applied (displayed in blue)

8.2.2 Resolution

Navigation: Home > Method Definition > Method name > Result and value handling > Resolution

Traditionally, moisture analyzers are operated at a maximum resolution of 1 mg / 0.01 %. The following setting allows you to determine moisture content using the resolution of the built-in analytical balance, which is ten times higher (0.1 mg weight / 0.001 % moisture). The high resolution is especially suitable for samples with low moisture content (lower than 1 %). Careful sample preparation is extremely important for these kind of samples!

Note

High resolution is not available for the display mode **ATRO moisture content** or **ATRO dry content**, **see** Settings of Display Mode (page 69).

You can define the following parameters:

Parameter	Explanation	Values
Resolution	Defines the readability of the built-in analytical balance.	Standard* High

* Factory setting

8.3 Workflow Handling

Navigation: Home > Method Definition > Method name > Workflow handling

With this functions you can define the workflow for the measurement

You can define the following parameters:

Menu item	Explanation	Further Information
Starting mode	Defines how the sample chamber is operated.	see Starting Mode (page 75)
Preheating	Defines the settings for preheating the sample	see Preheating (page 75)
	chamber before starting the measurement.	

8.3.1 Starting Mode

Navigation: Home > Method Definition > Method name > Workflow handling > Starting mode

In this menu item you can choose whether the sample chamber should operate automatically or manually (e.g. tare, stop drying).

You can define the following parameters:

Parameter	Explanation	Values
Starting mode	Defines how the sample chamber is operated.	Automatic*
		Manual

^{*} Factory setting

Automatic

Your instrument is set in the factory to the automatic operating mode. This mode can be used for most kind of samples. When closing the sample chamber, the weight of the sample is recorded and the measurement is started.

Manual

We advise you to use the manual operating mode for samples which contain readily volatile substances. In contrast to the automatic operating mode, in the manual operating mode the sample chamber does not automatically close when the [**Start drying**] button is tapped. However, the initial weight (wet weight) important for the determination of the moisture content is recorded. In the manual operating mode you have time for further preparation of the sample (e.g. mixing with quartz sand or even distribution of the sample) while weight losses due to evaporation during the preparation time are measured from the outset. As soon as the sample is ready for drying, press [‡] key. The automatic sample chamber closes and drying starts. In the manual operating mode you can open the sample chamber during a drying operation. In contrast to the automatic mode, drying will not be stopped but simply interrupted until the automatic sample chamber is closed again.

8.3.2 Preheating

Navigation: Home > Method Definition > Method name > Workflow handling > Preheating

If necessary, you can activate preheating function to preheat the sample chamber. Because of the short heat-up times with halogen moisture analyzers, no preheating is usually required for standard applications. A preheated instrument can improve the reproducibility and accuracy of the results, since the instrument is in the same climatic state with every measurement. Preheating will only executed until time of measurement (**Manual**), in standby mode even after the measurement.

You can define the following parameters:

Parameter	Explanation	Values
Preheating	Defines the settings for preheating the sample chamber (standby temperature) before starting the measurement.	OFF* ON Manual
	OFF = no preheating applied.	Standby

^{*} Factory setting

Manual

Before a measurement can be started, the instrument prompts the user to activate the function preheating. The instrument indicates, when it is enough preheated to start the measurement. The measurement can also be performed during the preheating phase by tapping **Stop preheating**. In this case, the duration of preheating will be recorded. After a maximal preheating time of 1 hour, preheat will be off.

You can define the following parameters:

Parameter	Explanation	Values
Temperature	Defines the standby temperature.	40100 °C (40 °C)*

^{*} Factory setting

Standby

The instrument indicates, when it is enough preheated to start the measurement. The measurement can also be performed during the preheating phase by tapping **Stop drying**. In this case, the duration of preheating will be recorded.

The function **Duration** allows you to deactivate preheating at a preselected time period after measurement. The feature **Turn-off time** allows to disable preheating at a defined clock time.

You can define the following parameters:

Parameter	Explanation	Values
Temperature	Defines the standby temperature.	40100 °C (40 °C)*
Duration	Defines the time which the instrument is maintained at standby temperature for a limited amount of time after measurement (time out).	Hours Minutes (3 h)*
Turn-off time	Defines the clock time at which Preheating will be off.	OFF* Hours : Minutes

^{*} Factory setting

8.4 General Method Properties

Navigation: Home > Method Definition > Method name > General method properties

8.4.1 Method name

Navigation: Home > Method Definition > Method name > General method properties > Method name You can define the following parameters:


Parameter	Explanation	Values
Method name	Renaming a method. Name should be unique and unambiguous.	any

8.5 Test a Method

Test a method

This function allows you to test the settings during the definition phase of a method. Testing is possible at any time. Test measurements are recorded in the journal and marked as test results. If you wish to work with the switch-off criterion «weight loss per unit of time» and are not familiar with the behavior of the sample, the test measurement will help you select the suitable setting. For more information about switch-off criteria, **see** Settings of Switch-off Criterion (page 67).

The diagram illustrates the progress of a drying. The points at which the individual switch-off criteria were reached (1-5) are marked.

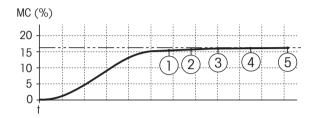

During this measurement, the measured values are recorded the switch-off criteria 1 to 5 and maybe one of a free-defined switch-off criterion. The measurement terminates after the defined time (default 30 minutes). Note that the start weight affects the switch-off criterion. The test time can also be switched off. In this case the measurement terminates after reaching the predefined switch-off criterion. Nevertheless, all switch-off criteria are recorded.

By comparing the measured values with values of a reference method e.g. drying oven and the standard deviation, you can define the appropriate parameter settings. For more information see the application brochure «Guide to Moisture Analysis».

Carry out all measurements under **Home > Method Definition > Method name > Tests** before releasing a method. The results of test measurements are specifically marked.

Attention

Before releasing the method, check if the determined switch-off criterion is set.

Performing a Test measurement

You perform a test measurement just like any other measurement. How to perform the measurement **see** Measurement (page 80).

Changing the test time

- Work screen for testing is displayed.
- 1 Tap **Test time** in the work screen.
 - ⇒ **Test time in min** appears.
- 2 Confirm with OK.

Displaying the parameter

- Work screen for testing is displayed.
- Tap the parameter in the work screen.
- ⇒ The parameter list appears.

Printing results of the determined switch-off criteria

For a printout example, See Information on Printouts (page 86) section «Special events».

Displaying results of the determined switch-off criteria

See Graphical Evaluations of Measurement Results (page 84)

9 Measurement

This function offers you to perform a measuring using previously defined methods. After choosing the measurement method, the measurement process can be started. The work screen guides you step by step through the measuring process.

The following functions are available:

- On the work screen, you can create a shortcut. This allows you to start a method directly from the home screen.
 - See Using Shortcuts (page 82).
- After completion of a measuring you can perform a further measuring with same method by tapping [Next Sample] or with an other method by tapping [Measurement].
- For graphical evaluations of measurement results tap [Results]. If you have defined the control limits in the method, passed, warned or failed status is shown.
 See Results (page 84).
- A detailed overview of the method parameter can you call up by tapping the parameter panel.
 See Work Screen (page 21).

9.1 Performing a Measurement

You are now familiar with all the parameters of your instrument and have defined all values for your sample. The instrument is now ready for the determination of your samples. In this Section you will learn how to perform measurements and how to stop the measurement process.

Switching on

- ► The instrument must be connected to the power supply for approximately 60 minutes in order to reach operating conditions.
- 1 To switch on the Instrument, press [也].
- 2 Log-in with your password, if needed.

Selecting method for measurement

- 1 Tap Measurement.
 - ⇒ The methods list is displayed.
- 2 Select the method to determine your sample.
 - ⇒ The work screen of the method appears.
 - ⇒ The sample chamber opens automatically.


Placing sample pan

- ► The display now prompts you to load the empty sample pan and tare the balance.
- 1 Place the empty sample pan in the sample pan handler.
- 2 Place the sample pan handler in the sample chamber. Ensure that the tongue of the pan handler fits exactly in the slot of the draft shield element. The sample pan must lie flat in the pan holder.

Note

We advise you to work with the sample pan handler at all times. The pan handler is ergonomic, automatic positioning, safe and provides protection against possible burns due to the hot sample pan.



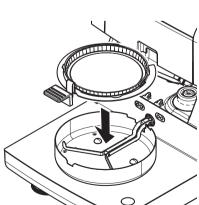


Taring balance

- Tap [->0/T<-].
 - ⇒ The sample chamber closes automatically for taring.
 - ⇒ When taring is completed, the sample chamber opens automatically.



Starting measurement

- After taring, the display prompts you to add the sample to the sample pan.
- 1 Add the sample to the sample pan. If you have defined a start weight, weigh the sample using the weighing-in aid.
- 2 Tap [Start drying].
 - ⇒ The sample chamber closes automatically.
 - ⇒ The drying process starts automatically.

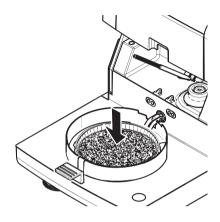

Drying process

You can follow the measurement process on the display, see Work Screen (page 21).

- The drying process is continuously displayed graphically.
- The current temperature in the heating module is displayed as well as the elapsed drying time and the current drying value.
- The display shows the selected settings.
- The drying process can be aborted by tapping [Stop drying].

At the end of the drying process you can read the moisture content of your sample on the display. If you have defined the control limits in the method, passed, warned or failed status is shown.

Removing Sample



CAUTION

Danger of burns


Sample, sample pan and sample pan holder may still be hot.

- ▶ The drying process is completed.
- ▶ The sample chamber is open (opens automatically).
- Carefully remove the sample pan handler from the sample chamber.

Note

To remove the sample pan from the handler, lift the pan slightly from below and remove it from the handler.

- To perform a further measurement with the current method, tap [Next Sample].
 - To perform a measurement with a new method, tap [Measurement].
 - To return to the home screen, press [1].

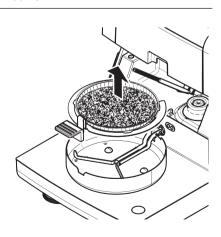

Stop drying

When you stop the measuring process by tapping [Stop drying], you can chose between two possibilities:

Abort without saving

Abort without saving of any data collected until now.

Abort measurement and save data

Already sampled data are stored and an entry will be made in the results. The result is marked as aborted.


Adding a comment

At the end of a measurement you can add a comment to the result of the measurement. This comment is stored in the measurement result and can be printed. A comment can only be entered before leaving the current measurement

- 1 To create a comment, tap [Note].
 - \Rightarrow The keyboard dialog appears.
- 2 Enter a comment.
- 3 Confirm with [**OK**].

9.2 Work with Weighing-in Aid

The weighing-in aid can be defined for each method and facilitates weighing in of the sample to a target value. This is particularly necessary if you require all the samples of a method you wish to process to have same weight in order to improve the repeatability of the measurement results. In addition, the weighing-in aid can be set as active so that the drying process cannot be started if the weight of sample is outside the set tolerance. You are therefore compelled to weigh-in the correct quantity of sample. If all sample weighed are within tolerances then this will improve the repeatability. The weighing-in aid is available only if start weight has been activated. For further information **see** Settings of Start Weight (page 71).

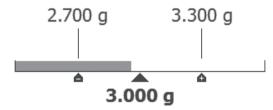

Icon	Function
	Lower weight limit (tolerance range)
	Target weight
A	Upper weight limit (tolerance range)

9.3 Using Shortcuts

Navigation for managing shortcuts: Home > Measurement > Method name > [7]

Shortcuts allow you to start methods directly from the home screen. Shortcuts are user-specific, i.e. each individual user can have his own shortcuts for the most common tasks.

You can define the following parameters:

Parameter	Explanation	Values
Shortcut name	Defines a name for the shortcut. It is recommended to choose a short name because only approx. 8 characters are displayed on the shortcut. Name should be unique and unambiguous.	any
Method name	Shows the name of the method. The name of the method is the none	
	target of the shortcut and can not be changed.	

Defining a shortcut

- Menu Measurement is activated.
- The method is selected.
- 1 Tap [7].
 - ⇒ Window **My shortcuts** appears.
- 2 Tap Add a shortcut to my home for this method.... (To abort, tap [X].)
 - ⇒ **New shortcut** appears.
- 3 To enter a name for the shortcut if needed, tap Shortcut name.

Note

It is recommended to choose a short name because only approx. 8 characters are displayed on the short-cut. Name should be unique and unambiguous.

- 4 Confirm with [**OK**].
- 5 To store the setting, tap [Save]. To abort, tap [Cancel].
- ⇒ The shortcut is added to the home screen.

Editing a shortcut

- ▶ The method is selected.
- 1 Tap [7].
 - ⇒ Window **My shortcuts** appears.
- 2 Tap Edit this shortcut.... (To abort tap [X].)
 - Shortcut properties appears.
- 3 To edit the name of the shortcut, tap **Shortcut name**.
- 4 Confirm with [OK].
- 5 To store the setting, tap [Save]. To abort, tap [Cancel].

Removing a shortcut


- ▶ The method is selected.
- 1 Tap [7].
 - ⇒ Window **My shortcuts** appears.
- 2 Tap **Edit this shortcut...** (To abort, tap [X].)
 - ⇒ **Shortcut properties** appears.
- 3 To remove the shortcut, tap [Remove].
 - ⇒ A message window appears.
- 4 Confirm with [Remove]. (To abort, tap [Cancel].)
- ⇒ The shortcut is removed from the home screen.

10 Results

10.1 Graphical Evaluations of Measurement Results

Navigation: Home > Results

This function allows you to manage and evaluate measurement results.

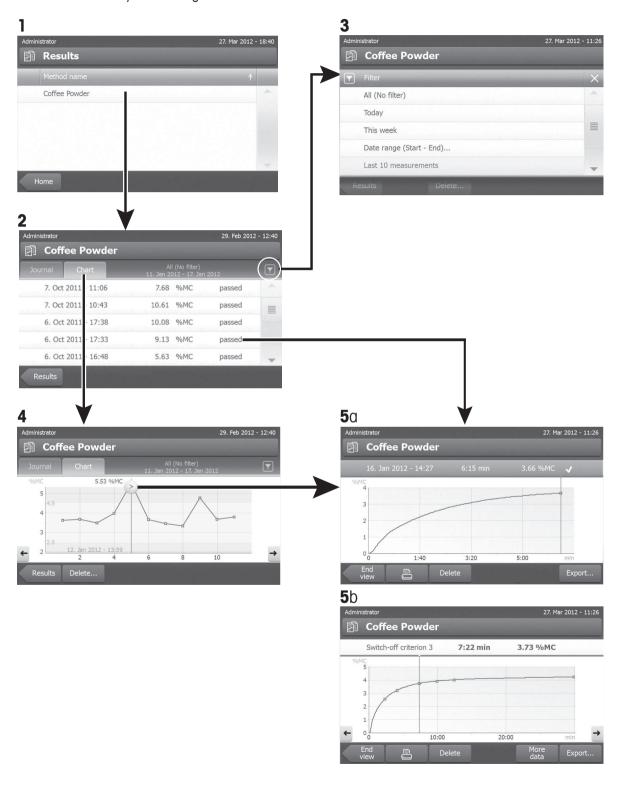

1 Method list

- Tap the method you want to evaluate.
 - ⇒ The journal view appears.

2 Journal view

The journal view allows you to start different graphical evaluations from a measurement series. You can perform the following functions:

- To start filter function, tap [\(\bigvarepsilon\)].
 - ⇒ The filter menu appears.
- To start chart view, tap [Chart].
 - ⇒ The chart view of the series of measurement appears (4).
- To start graphic view, tap a measurement result.
 - \Rightarrow The graphic view appears (5).

3 Filter menu

The filter function allows you to evaluate the measurement series according to various criteria. You can select the following criteria:

- All (No filter)
- Today
- This week
- Date range (Start End)...
- Last 10 measurements*
- Last 20 measurements
- Tap [X] to close filter menu.
- * Factory setting

4 Chart view

This view allows you to display the results of a series of measurements depending on the filter settings. If you have defined the control limits in the method, they are diagramed in the results.

The > marked measuring point is displayed with date, time, and the measurement result. You can perform following functions:

- To jump forward to the next measurement result, tap [->].
- To jump backwards to the previous measurement result, tap [<-]. Alternatively, tap directly on the desired
 measurement point.
- To recall the full measurement curve of the corresponding measurement, tap [>].

5a Graphics view

With this function you can display a graphical view of the detailed results of a single measurement. If you have defined the control limits in the method, passed, warned or failed status is shown, **see** Control limits (page 72).

You can perform the following functions:

Printing result

To print the result, tap [=].

Exporting results

See Exporting Results (page 87).

Deleting result

To delete this measurement result, tap [Delete] (depending on user rights).

More data

To display more measurement data, tap [More data]

5b Displaying results of test measurements

For test measurements **see** Test a Method (page 78).

This view allows you to display the results of the determined switch-off criteria. Each criterion is marked as a measuring point and displayed with its result. You can perform following functions:

- To jump forward to the next measurement result, tap [->].
- To jump backwards to the previous measurement result, tap [<-]. Alternatively, tap directly on the desired
 measurement point.

10.2 Information on Printouts

The illustrations in this chapter show examples of a standard measurement printout (factory setting) and a short measurement printout. The level of detail on the printouts depends on the selected settings in the menu.

Structure of the standard printout type

-MOISTURE DETERMINATION-METTLER TOLEDO Halogen MoistureAnalyzer HX204 SNR(Drying Unit) 2345 SNR(Terminal) 6788 SW(Drying Unit) 0.60.09 SW (Terminal) 4.20628 Method name TEST Drying prog Standard Drying temp 105°C Switch Off 3(1mg/50s) Display mode %MC Start weight OFF Control limits OFF Start mode Automatic User name Administrator Start weight 0.487 g 01:00 min 2.26 %MC 02:00 min 3.49 %MC Total time 0:22 min Dry Weight 0.470 g Moisture content 0.017 g End result 3.29 %MC Signature 14.10.2011 12:01 ----- END -----

Structure of the short printout type

on a character of the character type
-MOISTURE DETERMINATION-
METTLER TOLEDO Halogen MoistureAnalyzer
Type HX204 SNR(Drying Unit) 2345 SNR(Terminal) 6788 SW(Drying Unit) 0.60.09 Sw(Terminal) 4.20628
User name Administrator
Method name: TEST
Start weight 0.487 g 01:00 min 2.26 %MC 02:00 min 3.49 %MC Total time 0:22 min End result 1.79 %MC
14.10.2011 12:01
END

Special events

The **sample chamber was opened and closed** during drying process. When the sample chamber is open, drying is interrupted and then continued when the chamber is closed.

01:00	min	2.26 %MC
01:20	min	opened
01:28	min	closed
02:00	min	3.49 %MC

The **drying process has been aborted** by tapping [**Stop drying**] and the measurement result at the time of the abort is not recorded as it could be wrong.

01:00 02:00		2.26	
ABORTED			
14.10	.2011	12:01	L

Printing the test measurement

During the test measurement a record is printed which explains exactly when and with what measurement result each switch-off criterion was reached.

```
Switch-off crit. 3
Time 01:21 min
Dry weight 3.385 g
Moisture content
0.53 %MC
```

10.3 Exporting Results

Results can be exported to an external storage device e.g. memory stick. This data in CSV format can be imported into e.g. MS Excel for further evaluations.

Note

Results cannot be imported in a Moisture Analyzer.

You can define the following parameters:

Parameter	Explanation	Values
Location	Finds the location on the storage device.	Browse and select
File name	Defines the name of the new data file.	any
File type	Defines the type of the data file.	CSV*

^{*} Factory setting

Procedure

- Result is activated.
- ▶ An external storage device is connected e.g. memory stick.
- 1 To start, tap [Export]
 - ⇒ **Export Results** appears.
- 2 Tap Location > Browse and select.
 - ⇒ Location appears.
- 3 Select location for file and confirm with [OK].
- 4 Enter new file name, if needed.
- 5 To start export, tap [Export].

11 Maintenance



WARNING

Risk of electric shock

The instrument must be disconnected from the power supply, before cleaning or other maintenance work to be performed.

Note

- The thermal overload protection can not be reset by the user.
- The halogen lamp can not be replaced by the user.

In such cases, contact your METTLER TOLEDO representative.

11.1 Cleaning



CAUTION

Danger of burns

The interior parts of the heating module as well as the parts in the sample chamber could be very hot.

Wait until the heating module has cooled down completely.

To obtain precise measurement results, we recommend you to clean the temperature sensor and the protective glass of the halogen lamp regularly. Please note the following directions for cleaning your instrument.

General

Your Moisture Analyzer is made from high quality, resistant materials and can therefore be cleaned with a commercially available, mild cleaning agent e.g. isopropanol.

Attention

- Use a lint-free cloth for cleaning.
- Ensure that no liquids enters the interior of the instrument.

Heating Module

Clean the exterior of the heating module with a mild cleaning agent although the housing is extremely rugged and resistant to solvents.

• Terminal

On no account use cleaning agents, which contain solvents or abrasive ingredients, as this can result in damage to the terminal overlay.


 Never open the housing of the instrument – they contain no components, which can be cleaned, repaired or replaced by the user.

Note


After the temperature sensor or the protective glass have been cleaned, we recommend adjusting the heating module using the temperature adjustment kit, **see** Temperature Adjustment (page 54).

11.1.1 Sample Chamber

- ► Sample chamber is open.
- 1 Remove the sample pan handler (1), the sample pan holder (2) and the draft shield (3) for cleaning.
- 2 Carefully remove any deposits from the black temperature sensor (4).

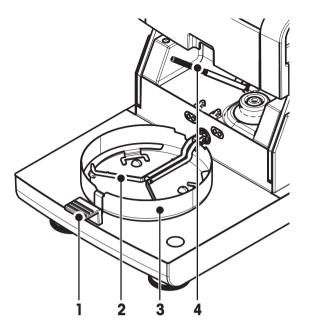

11.1.2 Heating Module

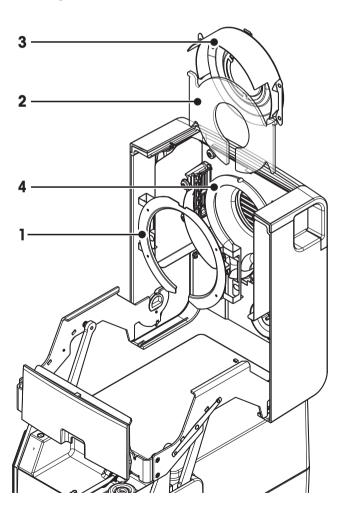

- 1 Reflector ring
- 2 Protective glass
- 3 Reflector with inspection window glass
- 4 Halogen lamp

To clean the protective glass, the reflector as well as the reflector ring, you should first open the heating module.

Attention

Avoid touching the round halogen lamp. If you have to remove any splashes, deposits, or spots of grease from the halogen lamp, it is advisable to use a weak organic solvent such as, for example ethanol. Ensure that the lamp is cool. **Do not remove the halogen lamp!**

Opening heating module for cleaning

- ▶ Sample chamber is open.
- 1 Inside, there is a locking device on each side. Push both outwards (together) for unlocking.

Note

Do not hold the module while unlocking it.

- ⇒ The upper part has been unlocked on both sides.
- 2 Open the heating module.




Removing reflector ring for cleaning

Note


For cleaning the protective glass, removing the reflector ring is not mandatory.

- Heating module is open.
- 1 Simultaneously, press the two lugs on both sides together.
- 2 Unhook the ring and remove it from the bracket.

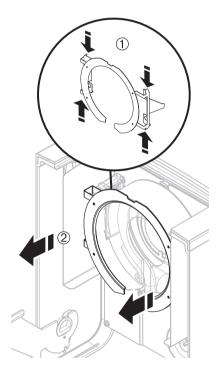


Removing reflector with inspection window glass for cleaning

- ▶ Reflector ring is removed.
- 1 To unlock, pull the spring clip forwards.
- 2 Pull out the reflector upwards off the bracket.

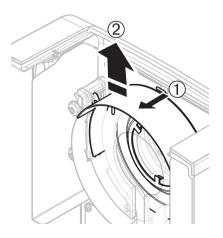

Removing protective glass for cleaning

- Reflector is removed.
- Pull out the protection glass upwards off the bracket.

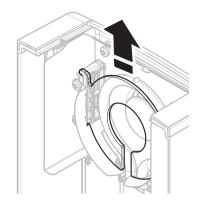

Reassembling after cleaning

Reassemble all parts in the opposite order.

- ► All parts are cleaned.
- 1 Insert the protective glass.
- 2 Insert the reflector with inspection window glass (until it clicks).
 Note


The lamp must be located in front of the reflector. Do not touch the lamp with your fingers.

3 Insert the reflector ring.

Attention

Observe the correct position! Assure that the ring is properly fitted.

4 Close the heating module (until it clicks).

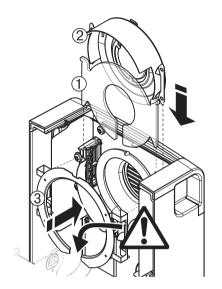


11.1.3 Fan Grill

The air inlet of the fan is located at the rear of the instrument and its exterior should be cleaned from time to time to free it from any dust deposits.

11.2 Dust Filter

If you use the Moisture Analyzer in a dusty environment with an dust filter, check the filter at regular intervals. Replacement Filters **see** Accessories and Spare Parts (page 100).

Replace the filter, if needed.

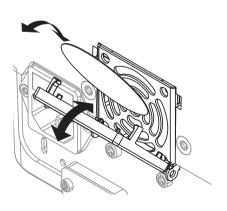

11.3 Replacing Power Line Fuse



CAUTION

Safety risk or damage of the instrument

Do not use a fuse of a different type or rated value, or short out (bridge) the fuse, because this can put your safety at risk and damage the instrument!

If the display of your terminal remains dark after switching on, in all probability the power line fuse of the drying unit is blown.

The power line fuse is located on the back of the drying unit. To change the fuse, proceed as follows:

- 1 Pull out the power cord.
- 2 Slide out the fuse holder (1) with a suitable tool such as a screwdriver.
- 3 Remove the fuse (3) and check its condition.
- 4 If the fuse is blown, replace the fuse with one of the same type and the same rated value (5 x 20 mm, T6.3H 250 V).

Note:

A spare fuse is located in the holder (2).

Fuse **see** General Technical Data (page 97)





In conformance with the European Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

Thank you for your contribution to environmental protection.



12 Troubleshooting

Errors can occur during operation of your instrument. This section describes how you can rectify these errors.

12.1 Error Messages

Most error messages appear in plain text directly in the respective application, and usually accompanied by a text describing how to correct the error. Error messages of this type are self-explanatory and therefore are not mentioned below. The following error messages can appear instead of the weighing result.

Error Message	Cause	Rectification				
Weight display						
	Overload — The weight on the pan exceeds the weighing capacity of the balance.	Reduce the weight of the sample.				
	Underload – Sample pan holder is missing.	Insert sample pan holder. If needed, reboot the system by disconnecting and connecting to the mains.				
0.0000	Weight display flashes / Out of zero range — When the instrument was switched on or upon zeroing, one or more limits were exceeded. The usual reason for this message to appear is when there is a weight on the weighing pan when the balance is switched on.	 Remove the weight. 				
Adjustment	I					
Unstable weight.	No stability during adjustment.	 Ensure ambient conditions and an optimum location. Take care that no part of the sample or the sample pan touches the draft shield or the sample pan handler. Ensure that the sample pan holder is correctly installed as well as not defect. Highly volatile substances in the sample also prevent a stable weighing result being detected since the sample is continuously losing weight. 				
Weight out of range.	Either no weight , or the wrong weight , has been placed on the sample pan during adjustment. (This message is also displayed if you do not remove the weight when prompted to do so by the instrument.)	Repeat the adjustment process and load the required adjustment weight.				

12.2 Status Messages

Status messages are displayed by means of small icons in the status bar of the display. For further information **see** Status Message Icons (page 18). The status icons indicate the following:

	Cause	Rectification
<u>\$\$\$</u>	Hot surface Indicates that the temperature inside the sample chamber is over approximately 50°C. The parts in the sample chamber and the sample itself can be very hot and there is a risk of burns.	The status icon will disappear as soon the temperature inside the sample chamber is less than approximately 50°C.
©	The built-in level sensor has determined that the instrument is not correctly leveled.	 Level the instrument immediately. See Leveling Drying Unit (page 26)
		⇒ If the instrument is leveled correctly, the icon disappears.
€	The instrument would like to carry out a fully automatic FACT adjustment but is unable because another operating sequence is running.	The adjustment is carried out as soon as the balance is unloaded, the display becomes stable and no key has been pressed for 2 minutes. The status icon disappears after the adjustment is successfully completed.
<u></u>	The battery in your instrument must be replaced. This battery ensures that the date and time are not lost when the instrument is disconnected from the power supply.	 Contact your METTLER TOLEDO representative as soon as possible. ⇒ A service technician will change the battery.
Se	Your instrument is due for servicing.	Contact your METTLER TOLEDO representative as soon as possible to have a technician service your instrument.

12.3 What to Do if...

Symptom	Countermeasure
Display remains dark after switching on	 Ensure that the terminal is properly connected to the drying unit. Ensure that the instrument is connected to the power supply and the power is on. Check the power line fuse of the drying unit and replace if necessary, see Replacing Power Line Fuse (page 92). If the problem persists, contact a METTLER TOLEDO representative.
Keys and buttons do not respond	 Reboot the system by disconnecting and connecting to the mains. If the problem persists, contact a METTLER TOLEDO representative.
Printer that is connected does not print	 Ensure that the printer is on and enabled in the menu. See Peripherals (page 34) Check the printer settings. See Recommended Printer Settings (page 105).
Incorrect characters are printed	 Change the bit/parity setting of the printer and the instrument to 8/NO. Check if both instruments have the same baud rate setting, see Peripherals (page 34). Use the correct character sets, see Recommended Printer Settings (page 105).
Measurement takes too long	 You have selected an unsuitable switch-off criterion, see Settings of Switch-off Criterion (page 67). An excessive amount of sample can also be the cause of slow drying, likewise samples which tend to form a skin which hinders vaporization. Perform a measurement at higher temperature. Enlarge the surface of the sample, e.g. by crushing or grinding. Use absorbent glass fiber filters for liquids. If the sample is very temperature sensitive and decomposes, reduce the temperature. If measuring is unstable, check correct positioning of pan, sample handler, sample, draft shield or sample holder.
Instrument does not heat after the start	Halogen lamp is defective or the heating module is overheated and the thermal overload protection has switched off the heating. In case, contact your METTLER TOLEDO representative.

Symptom	Countermeasure
Measurement results are not repeatable	Support on which the instrument is standing is not sufficiently stable. Use a stable support.
	• Surroundings are very unstable (e.g. vibrations, air draft, humidity). Provide better environmental conditions.
	Sample has more or less moisture between sampling and starting the drying process.
	Sample is not evenly dispersed on the pan.
	Start weight has not always the same value.
	 Samples are not homogeneous, i.e. they have different compositions. The more inhomogeneous a sample, the larger the amount of sample needed to obtain a repeatable result.
	 Selected drying time is too short for the "Timed switch-off" criterion. Extend the drying time or select a suitable switch-off criterion "Weight loss per unit of time".
	• Sample does not become completely dry (e.g. awing to skin formation). Dry the sample with the aid of glass fiber discs.
	Selected temperature is too high and the sample has oxidized or decomposed. Lower the drying temperature.
	Sample boils and the splashed drops continuously change the weight. Lower the drying temperature.
	Granulation is not homogeneous or too large.
	Insufficient heating power because the protective glass of the halogen
	lamp is dirty. Clean the protective glass, see Cleaning (page 88).
	Temperature sensor is contaminated or faulty. Clean the temperature sensor, see Cleaning (page 88).
	If the problem persists, contact a METTLER TOLEDO representative.

13 Technical Data

13.1 General Technical Data

Drying Unit

Heating Module Halogen ring-shaped radiator

Temperature range 40–230 °C

Temperature step 1 °C

Temperature programs standard, rapid, gentle, step

Balance

Maximum load 200 g

Readability 1 mg / 0.1 mg

Minimum sample weight 0.1 g
Weighing technology Monobloc

Adjustment FACT, internal weight, external weight

Moisture Content

Readability 0.01%
Readability, fine range 0.001%
Repeatability (sd) with 2 g sample max. 0.05%
Repeatability (sd) with 10 g sample max. 0.01%

Materials

Drying Unit

Housing Plastic, PBT, Crastin S0653-GB20 Inspection window grill Plastic, PEEK-HT G22 (UL94-V0)

Protective glass Glass ceramics Halogen lamp Quartz glass

Reflector Stainless steel, X2CrNiMo17-2 (1.4404)

Reflector bracket Plastic, PEEK-HT G22 (UL94-V0)

Draft shield, interior bottom plate Stainless steel, X2CrNiMo17-2 (1.4404)

Terminal

Top housing EN ZL-ZnAl4Cu1 (EN ZI-0410)

Bottom housing PA12 GB30

Protection and Standards

Overvoltage category Class II
Degree of pollution 2

Standards for safety and EMC see Declaration of Conformity (part of standard equipment)

Range of application for use in dry interior rooms

Environmental Conditions

Height above sea level up to 4000 m Ambient temperature range 5 °C to 40 °C

Relative air humidity 10% to 80% at 31 °C, linearly decreasing to 50 % at 40 °C,

noncondensing

Warm-up time At least 60 minutes after connecting the instrument to the power

supply; when switched on from standby, the instrument is ready

for operation immediately.

Power Supply

110 V AC Version 100 V–120 V, 50/60 Hz, 4 A 230 V AC Version 220 V–240 V, 50/60 Hz, 2 A

Voltage fluctuations -15%+10%

Power load max. 450 W during drying process

Power line fuse 5 x 20 mm, T6.3H 250 V

Interfaces

Drying Unit

1x System (Terminal – drying unit)

1 x RS232C (9-pin socket)

2 2 V V P V Last (Torre A socket)

2x USB Host (Type A socket)

1x Memory Card slot SD/SDHC up to 32 GB (SDXC not sup-

ported)

Data

Export data format of results *.CSV

Comma-separated values (CSV) file

Hardware

Drying unit

Sample chamber opening/closing Motorized

Leveling 2 leveling screws,

Level indicator and Inclinometer

Sample pan Ø 90 mm Maximum heigh of sample 15 mm

Thermal overload protection Bimetallic-element switch in heating module

Dimensions (w x h x d) 199 x 139 x 428 mm

See Dimensions (page 99)

Weight, ready to measure 6.8 kg

Terminal

Display WVGA color display, touch-sensitive

Reading angle Adjustable, 2 folding feet

Dimensions (w x h x d) 200 x 63.5/79.5 x 134.5 mm


See Dimensions (page 99)

Weight 1.2 kg


13.2 Dimensions

(all dimensions in mm)

Drying Unit

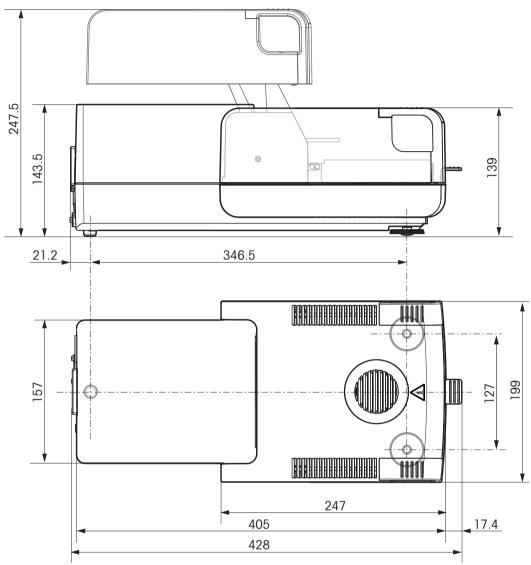

Terminal

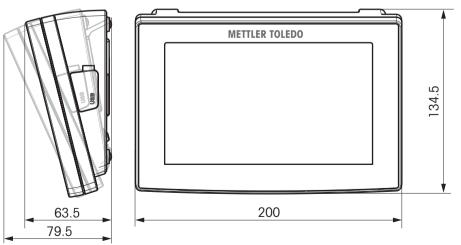

14 Accessories and Spare Parts

Description Part No.

Power supplies



Country-specific 3-Pin power cable with grounding conductor.

Power cable CH	00087920
Power cable EU	00087925
Power cable US	00088668
Power cable IT	00087457
Power cable DK	00087452
Power cable GB	00089405
Power cable AU	00088751
Power cable ZA	00089728
Power cable BR	30015268
Power cable JL	00225297
Power cable IN	11600569
Power cable JP	11107881
Power cable TH, PE	11107880

Cables for RS232C Interface



RS9 - RS9 (m/f): connection cable for PC, length = 1 m

11101051

Cables for USB Interface



USB (A -B) connection cable for connection to PC, length = 1 m

12130716

Cables for Terminal



Terminal cable, length = 0.68 m


30003971

Printers



RS-P25 printer with RS232C connection to instrument	11124300
Paper roll, set of 5 pcs	00072456
Paper roll, self-adhesive, set of 3 pcs	11600388
Ribbon cartridge, black, set of 2 pcs	00065975

	RS-P42 printer with RS232C connection to instrument Paper roll, set of 5 pcs Paper roll, self-adhesive, set of 3 pcs Ribbon cartridge, black, set of 2 pcs	00229265 00072456 11600388 00065975
Sample Pans		
	Aluminum sample pan, HA-D90, set of 80 pcs	00013865
	Professional aluminum sample pan, extra strong, set of 80 pcs	11113863
	Stainless steel reusable sample pan 6 mm, DA-DR1, set of 3 pcs	00214462
Adjustment Parts	Certified adjustment weight, 100 g (F1)	11119531
	OIML / ASTM Weights (with calibration certificate) see www.mt.com/weights	ghts
	HX/HS Temperature adjustment kit, certified	30020851
	SmartCal™, Moisture Analyzer Reference Substance	
The state of the s	cSmartCal™, certified, 12 tests	30005793

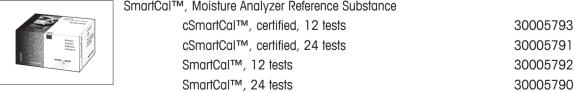


Anti-theft Devices



Steel cable

11600361

Miscellaneous

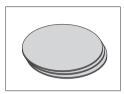


Glass fiber filter (for Liquids), set of 100 pcs

00214464



Dust filter, set of 50 pcs

30020838

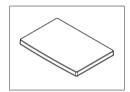


Protective cover for Terminal

30003957



Transport case


30020836

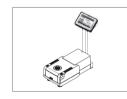

Stand for terminal


30018474

Spare Parts

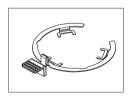


Sample pan handler

30020852

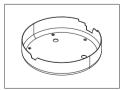

Draft shield

30007150

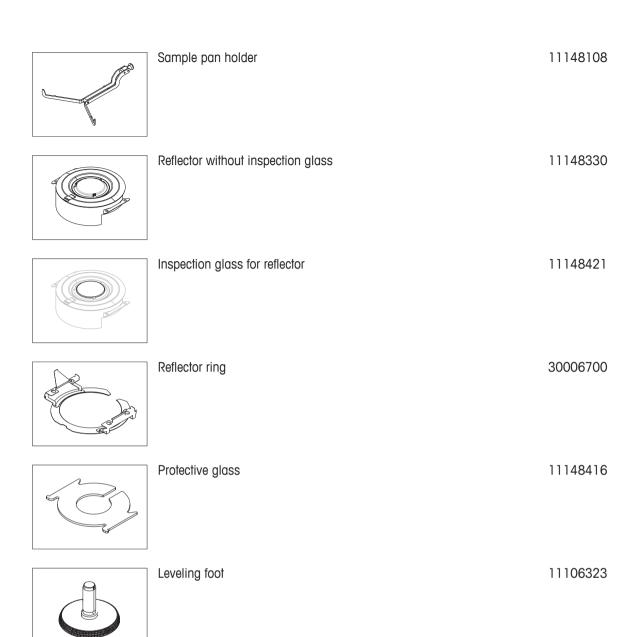


15 Appendix

15.1 How to Obtain Best Results

In this Section you will find important information on how to obtain optimum results. You will discover which parameters influence the measurement process and how you can optimize the settings of your instrument to achieve the best measurement results.

15.1.1 Measurement Principle of Halogen Moisture Analyzer

Your instrument performs measurements based on the **thermogravimetric principle**, i.e. the moisture is determined from the weight loss of a sample dried by heating.

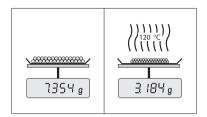

In principle, your instrument thus comprises two instruments: An analytical balance and a heating module. In contrast to other thermogravimetric methods (drying oven, infrared, microwave), the Halogen Moisture Analyzer operates with halogen heating technology. This ensures fast heating of the sample and thus guarantees rapid availability of the measurement results.

Irrespective of the measurement method, a correct preparation of the sample and the correct choice of the following measurement parameters improve the quality of the measurement result:

- Sample size
- Drying temperature
- Switch-off mode
- Drying time

Attention

Inappropriate setting of these parameters can cause the results to be incorrect or misleading. For this reason, check that the results for each type of sample are what might reasonably be expected.

You will find detailed information on the relationships between these parameters in the delivered application brochure «Guide to Moisture Analysis» or **see** Accessories and Spare Parts (page 100).

In practice not only the quality of the measurement results, but also the speed of the measurement process is important. Thanks to its drying principle (heat generated by a halogen radiator), the Halogen Moisture Analyzer is very fast. You can increase the speed even further through optimum setting of the instrument, e.g. using the **Drying programRapid**.

The optimum drying temperature and the drying time are dependent on the nature and size of the sample and on the desired accuracy of the measurement results. These can be determined only by experiment. The Halogen Moisture Analyzer supports you in this task: It offers result recording of test measurements in the menu **Method Definition**.

15.1.2 Notes on Adjustment of Balance and Heating Module

The balance and the heating module in your drying unit can be adjusted using the appropriate accessories, **see** Accessories and Spare Parts (page 100). The heating module and the balance can also be tested (to check the adjustment). The user can define a test weight or the test temperature, together with permissible tolerances for this test. A test report can be printed out showing the test result has passed or failed.

A Moisture Analyzer is typically used in place of or in addition to the drying oven method. In an oven heat energy is transferred by the flow of air, which establishes an equilibrium between the sample temperature and the ambient temperature. This is not the case in a Moisture Analyzer. The actual sample temperature primarily depends on the specific absorption properties of the sample (dark samples absorb more heat), which can change during the measurement process. There can also be differences between the temperature at the surface of the sample and the temperature inside the sample. The heat output is therefore not dependent on the true sample temperature but instead is regulated by a temperature sensor underneath the halogen heating module.

For the reasons explained above, the sample temperature will be slightly different than the temperature shown on the instrument display. By regularly testing or adjusting the drying unit, you will ensure a consistent and reproducible heat output for the entire lifetime of your instrument.

Note:

- METTLER TOLEDO offers an adjustment service please contact your METTLER TOLEDO representative.
- We recommend to adjust the instrument exclusively under operating conditions.
- After the temperature sensor or the protective glass have been cleaned, we recommend adjusting the heating module using the temperature adjustment kit, see Accessories and Spare Parts (page 100). For the procedure for testing or adjusting the drying unit (balance/heating module), see Test/Adjust (page 52).

15.1.3 Optimum Sample Preparation

Preparation of the sample is decisive for the speed of the measurement process and the quality of the measurement results.

Note:

Basic rules for the preparation of your sample:

The amount of sample you select should be as small as possible and only as large as necessary.

Excessive amounts of sample require more time for drying and thus prolong the measurement process. If the amount of sample is too small, the measurement result may possibly not be representative of the true moisture content. The following always holds: The greater the inhomogeneity of the sample, the larger the amount of sample needed to obtain a repeatable result.

Distribute the sample evenly over the sample pan

You thus increase the surface area of the sample and facilitate heat absorption. The base of the pan should be evenly covered.

With liquid, fat-containing, melting and highly reflecting samples, you should use the sample with the glass fiber filter available as optional equipment, **see** Accessories and Spare Parts (page 100). This also applies to samples which form a skin on their surface when heated. The glass fiber filter ensures even and rapid heat distribution and prevents the formation of a skin on the sample surface.

15.1.4 Further Information on Moisture Determination

You will find further information on the moisture determination, the importance of the parameters and the preparation of the samples in the delivered application brochure «Guide to Moisture Analysis» dealing with moisture determination, **see** Accessories and Spare Parts (page 100).

Useful tips and a variety of example methods (comparison between Halogen Moisture Analyzer results and the oven drying method) can be downloaded from:

www.mt.com/moisturewww.mt.com/moisture-methodswww.moisture-guide.com

If you require information on specific applications, your METTLER TOLEDO Customer Service will be pleased to help you.

Attention

Moisture determination applications must be optimized and validated by the user according to local regulations. Application-specific data provided by METTLER TOLEDO is intended for guidance only.

15.2 Recommended Printer Settings

English, German, French, Spanish, Italian, Polish, Czech, Hungarian

Printer		•	
Model	Char Set	Auto baudrate	Balance feature
RS-P25	IBM/DOS	Off	Disable

Printer						
RS-P42	IBM/DOS 1)		-	_		_
Instrument / Printer						
Model	Baudrate	Bi	t / Parity	Stop Bits	;	Handshake
RS-P25	9600		8/NO	1		Xon/Xoff

8/N0

Xon/Xoff

Brazil Portuguese

RS-P42

Printer			
Model	Char Set	Auto baudrate	Balance feature
RS-P25	IBM/DOS	Off	Disable
RS-P42	2)	_	_

Instrument / Printer						
Model	Baudrate	Bit / Parity	Stop Bits	Handshake		
RS-P25	9600	8/NO	1	Xon/Xoff		
RS-P42	2)	2)	2)	2)		

¹⁾ Printer settings not available.

15.3 Third Party License/Notice

This section contains Third Party Software Notices and/or Additional Terms and Conditions for licensed third party software components included within SOFTWARE PRODUCT.

This SOFTWARE PRODUCT is based in part on the work of:

Qwt project

For user's guide **see** http://qwt.sf.net

For LGPL license V2.1 see http://www.gnu.org/licenses/old-licenses/lgpl-2.1.html

KompexSQLite

For LGPL license V3 see http://www.gnu.org/licenses/lgpl.txt

1200

• Qt library 4.8

For LGPL license V2.1 **see** http://qt-project.org/doc/qt-4.7/lgpl.html For GPL license V3 **see** http://gt-project.org/doc/qt-4.7/lgpl.html

decNumber

For ICU license V3.68 see http://source.icu-project.org/repos/icu/icu/trunk/license.html

• Simpleini

For MIT license **see** http://www.opensource.org/licenses/mit-license.php

²⁾ Required font for this language not available.

10	6 Index			Disposal Draft shield element	92 12
A	Access rights Accessories Account Policies User Adding comments Adjustment Balance FACT	44, 46 100 47 46 82 17, 28 52 52		Drying Steps Temperature Unit Unit dimensions Drying program Gentle Rapid Standard Dust filter	67 64, 65, 65 12, 25, 97, 98 99 65 64 63 12, 91
_	Heating module Notes Printouts Settings Touch screen Weight Weight external Weight internal Anti-theft Appendix ATRO Automatic sample chamber	52 104 56 40 36 53 53 53 12, 27 104 73	E	Eco Edit Method Shortcut Entering Numerical values Text and numbers Error messages Export Result Results Settings and methods	76 62 83 19 19 93 87 98 48
B C	Back up Basic elements Battery Brightness Chart	49 18 94 33, 36	F	FACT Fan Filter Firmware update First measurement Fuse	28, 52, 94 12, 91, 91 84 50 29 92, 98
	Chart view Cleaning Color Comments to measurement Commissioning Connecting the Instrument Connectivity Control limits Copy Method CSV format	22 88 33, 36 82 23 24 12, 14 72 62 87, 98	G	General method properties General safety information General settings Gentle drying Graphics panel Graphics view Groups	77 8 17 65 22 22, 84 44
D	Data Data management Date Default settings Define	98 48 20, 27, 34 35 62 83 62 99 33, 35 69	Ī	Halogen lamp Heating module History Home screen Host settings Identifications Identifications panel Import Settings and methods Inspection window Installation Intended Use	12 12, 12, 88 60 15, 17 34 38 22 48 12 23 8

	Interface	98	_	Overload	93
	Interim results	38	Р		
	Introduction	7		Parameter panel	22
J				Password	33, 46
	Journal	84		Peripheral settings	34
K				Power line fuse	12, 92
•	Kensington slot	12, 27		Power supply socket	12, 98
	Keyboard	12, 27		Power supply socket Preheating	12 75
	text and numbers	19		Printer	73
	Keyboard language	33		Settings	34, 105
	Keypad			Printout	,
	numerical values	19		Intermediate results	38
	Keys on the terminal	15		Language	38
L				Short Standard	86
	Languages	33, 35		Protective clothing	86 9
	Printout	38	_	1 Tolective clothing	
	Level		R		
	Indicator	12		Ramp time	66
	Sensor Leveling	12, 27, 94 26		Rapid drying	64
	Leveling screw	12		Reading angle	27
	Limits	72		Remove Method	62
	Lists	21		Shortcut	83
	Location	23		Replace	00
	Login	47		Dust filter	91
М				Power line fuse	92
•••	Maintenance			Reset	50
	Cleaning	88		Resolution Restore	73 49
	Dust filter	91		Result	49
	Heating module	88		Export	87
	Power line fuse	92		Result and value handling	72
	Measurement	17		Results	17, 84, 84
	First	29		Results export	98
	Measurement principle	104		RS232C	13, 14
	Memory Card Method	14	S		
	Сору	62		Safety Information	
	Definition	61		General	8
	Edit	17, 62		Intended use	8
	Export and import	48		Protective clothing	9
	Name	22, 77, 77		Signal words	8
	New	17, 62		Staff safety	9
	Properties Remove	77 62		Warning Symbols	8 12, 15
	Renaming	77		Sample chamber Sample pan handler	12, 13
	Test	78		Sample pan holder	12
	MT-SICS	34		Sample preparation	105
N				Scope of delivery	23
IN	AL II I	22		Screen settings	33, 35, 36
	New method	62		SD card	14
	Numbers Numerical values	19 19		Selecting the location	23
_	TAUTHELICAL VALUES	10		Setting up drying unit	25
0				Settings Adjustment	17 40
	Out of zero range	93		Aujusiiileiii	40

	Date	20, 27, 34		Temperature	58 79
	Export and import Host	48 34		Time Weight external	78 57
	Peripheral	34		Weight internal	57 57
	Regional	34		Text	19
	Screen	33			104
		34		Thermogravimetric Time	20, 27, 34
	System Temperature Test	41		Touch screen	20, 27, 34 15
	Temperature Test Time	20, 27, 34			36
	Weight adjustment	20, 27, 34 42		Touch screen adjustment Troubleshooting	93
	0 ,	41		Troubleshooling	
	Weight test Shortcuts	17, 18, 22, 82	U		
	Define	83		Underload	93
	Edit	83		Unpacking	23
	Remove	83		Update	51
	Show results	17		USB	13
	Signal words	8		USB Host	14
	Software update	50		User	
	Spare parts	100		Account	46
	Staff Safety	9		Groups	44
	Stand for terminal	26		Home screen	17
	Standard drying	63		Interface	16
	Standard equipment	23		Management	44
	Standby	15, 76		Profile see User manage-	44
	Start measurement	17		ment	
	Start weight	71, 82		Shortcuts	17, 82
	Starting mode	75	_		
	Startup user	47	V		
	Status icon	94		Value panel	22
	Status messages	18, 93	W		
	Stop drying	82, 87	•••	Warning Symbols	8
	Switch button	20		Weighing-in aid	71, 82
	Switch ON/OFF	15		Weight	/1,02
	Switch-off criterion	67, 78, 87		Adjustment external	53
	Switch-off time	76		Adjustment internal	53
	System	12, 14		Test external	57
	System management	48		Test internal	57 57
	System settings	34		Weight adjustment	0,
Т				Settings	42
•	Tables	21		Weight test	
	Technical data	97		Settings	41
	Temperatur	37		Work screen	21
	Adjustment settings	42		Workflow handling	75
	Temperature	42		<u> </u>	
	Adjustment kit	12			
	Sensor	12			
	Standby	76			
	Switch-off criterion	67			
	Test	58			
	Test settings	41			
	Terminal	27, 88			
	Terminal dimensions	99			
	Test	17			
	Balance	57			
	Heating device	57			
	Measurement	78, 87			
	Method	78, 87			

GWP® – Good Weighing Practice™

The global weighing guideline GWP® reduces risks associated with your weighing processes and helps to

- choose the appropriate balance
- reduce costs by optimizing testing procedures
- comply with the most common regulatory requirements

www.mt.com/GWP

www.mt.com/moisture

For more information

Mettler-Toledo AG, Laboratory & Weighing Technologies

CH-8606 Greifensee, Switzerland Tel. +41 (0)44 944 22 11 Fax +41 (0)44 944 30 60 www.mt.com

Subject to technical changes. © Mettler-Toledo AG 04/2012 30019565 en

