
<b>English</b>	<b>Operating Instructions</b> BT-P42 Printer
<b>Deutsch</b>	<b>Bedienungsanleitung</b> Drucker BT-P42
<b>Français</b>	<b>Mode d'emploi</b> Imprimante BT-P42
<b>Español</b>	<b>Instrucciones de manejo</b> Impresora BT-P42
<b>Italiano</b>	<b>Istruzioni per l'uso</b> Stampante BT-P42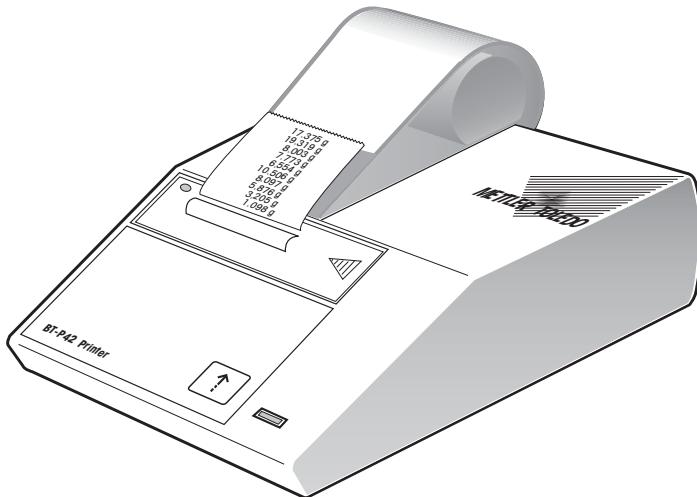

 **Bluetooth®** certified

Bluetooth is a registered trademark of Bluetooth SIG, Inc.

English

Deutsch

Français

Español

Italiano

**METTLER** **TOLEDO**

# Content

<b>1</b>	<b>Introduction .....</b>	<b>4</b>
<b>2</b>	<b>Startup Procedure .....</b>	<b>4</b>
2.1	Safety measures.....	4
2.2	Power supply.....	4
2.3	Inserting paper.....	5
2.4	Connecting the printer to the device .....	5
<b>3</b>	<b>Operation.....</b>	<b>5</b>
<b>4</b>	<b>Printer Test .....</b>	<b>6</b>
<b>5</b>	<b>Maintenance .....</b>	<b>6</b>
5.1	Changing paper .....	6
5.2	Replacing the ribbon.....	6
5.3	Cleaning.....	7
5.4	Disposal.....	7
<b>6</b>	<b>Malfunctions .....</b>	<b>7</b>
<b>7</b>	<b>Consumables and Spare Parts .....</b>	<b>8</b>
<b>8</b>	<b>Technical Data .....</b>	<b>9</b>

# 1 Introduction

## What are the capabilities of the BT-P42 Printer

The BT-P42 is a rugged and simple to operate dot matrix printer for METTLER TOLEDO balances with an a wireless Bluetooth® interface. The printout on normal paper fulfills the requirements of modern quality assurance systems (GLP, GMP, ISO 9001, etc.). These define, among other things, the principles and duties for the user of measuring instruments or inspection, measuring and test equipment. The BT-P42 complies with the principle of traceability of measurements by precisely documenting the origin of data, as well as calibration procedures. It fulfills the obligation to archive measurement results by its thermally stable and lightfast print. On connection to METTLER TOLEDO balances, the BT-P42 prints all data sent by the balance without restriction, that is not only weight values and calibration procedures, but also the results of all applications, built into the balance.

# 2 Startup Procedure

## 2.1 Safety measures

Always operate and use your product only in accordance with the instructions contained in this manual. The instructions for setting up your new product must be strictly observed. The BT-P42 has been tested for the connections and intended purposes documented in these Operating Instructions. However, this does not absolve you from the responsibility of performing your own tests of the product supplied by us regarding its suitability for the methods and purposes you intend to use it for. You should therefore observe the following safety measures.

- The BT-P42 Printer may be used only indoors and on no account in hazardous areas.
- The electronics of the BT-P42Printer are protected against the ingress of splashing water and dust. However, as the printer housing is not tight the printer should not be used in the presence of liquids.
- Comply with cleaning instructions (see item 5.3), protect power plug against wet conditions.
- Never open the printer housing. Should you ever have problems with your printer, please contact your responsible METTLER TOLEDO dealer.

## 2.2 Power supply

Your printer is supplied with an AC adapter with a country-specific power cable. The power supply is suitable for all line voltages in the range: 100 - 240 VAC, 50 - 60 Hz (for exact specifications, see item 7).

### Attention


- First, check the local line voltage is in the range 100 - 240 VAC, 50 - 60 Hz and whether the power plug fits your local power supply connection. **If this is not the case, on no account connect the printer or the AC adapter to the power supply,** but contact the responsible METTLER TOLEDO dealer.
- **Only plug the adapter into a socket which is grounded.**

### Important

- Before operating, check all cables for damage.
- Guide the cables so that they cannot become damaged or interfere with the weighing process!
- Take care that the AC adapter cannot come into contact with liquids!
- The power plug must be always accessible.


## 2.3 Inserting paper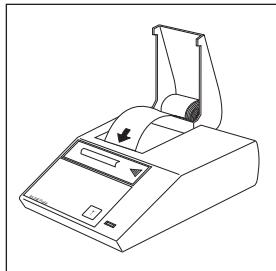

- Feed paper through the slot in the printing unit in the direction of the arrow.
- Switch on printer with toggle switch at the rear.
- Press and hold until sufficient paper has been fed in.

English

## 2.4 Connecting the printer to the device

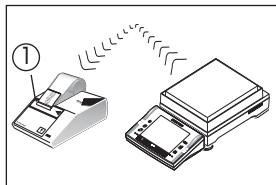- Position the printer within range of the balance.  
Maximum distance: 10 m
- Connect the device in accordance with the Installation and Operating instructions for the BT or BTS interface option.  
  
As soon as the device to be connected has been set up for the BT-P42, the printer is ready to operate. The connection may take 10 to 20 seconds to establish.  
  
The connection status is indicated by the blue Bluetooth lamp (1):
  - Lamp flashing: No connection
  - Lamp continuously lit up: Connection present

## 3 Operation

- Pressing the printer key advances the paper 1 line (line feed).
- Pressing the balance key with the symbol prints out the stable weight value.  
As long as the balance does not show a stable weight value, nothing will be printed out.

## 4 Printer Test

The BT-P42 is equipped with a self-test which automatically prints out the character set of the printer.

### Starting the test

- Switch on BT-P42 and at the same time press and hold for a few seconds.

### Terminating the test

- Switch off the BT-P42.


## 5 Maintenance

Under normal ambient conditions, the BT-P42 Printer requires practically no maintenance thanks to its rugged construction. The METTLER TOLEDO service department is available should servicing be required. Please ask your responsible METTLER TOLEDO dealer for details.


### 5.1 Changing paper

Insertion of a new roll of paper is described at item 2.4. Order number for paper, see item 7.

### 5.2 Replacing the ribbon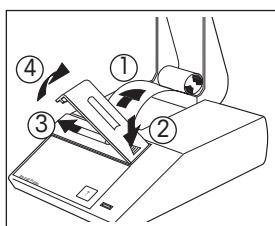


- Pull paper out of printing unit (1).
- Remove ribbon cover (2, 3, 4).
- To remove ribbon cartridge, press in direction of arrow (5).
- Insert new ribbon and if need be, take up slack with tensioning wheel (6).
- Refit ribbon cover.
- Insert paper and feed through slit in the printing unit as far as it will go.
- Press and hold until sufficient paper has been fed in.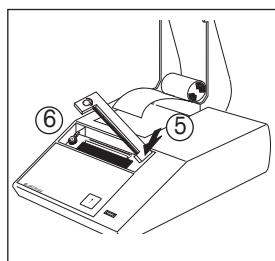

## 5.3 Cleaning

As the printer housing is made of top grade, resistant materials, all commercially available cleaning agents may be used.

## 5.4 Disposal

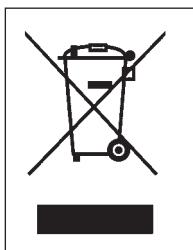In conformance with the European Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

Thank you for your contribution to environmental protection.

## 6 Malfunctions

Malfunction / Message	Possible cause	Rectification
Printout unreadable	<ul style="list-style-type: none"><li>Ribbon worn / entangled</li><li>Service life of printing unit at an end</li></ul>	<ul style="list-style-type: none"><li>Change / tension ribbon</li><li>Please contact METTLER TOLEDO service</li></ul>
Green pilot lamp does not light up	<ul style="list-style-type: none"><li>No power</li><li>fuse defective</li></ul>	<ul style="list-style-type: none"><li>Printer on</li><li>Please contact METTLER TOLEDO service</li></ul>
Blue Bluetooth display does not light up	<ul style="list-style-type: none"><li>No power</li><li>Bluetooth module faulty</li></ul>	<ul style="list-style-type: none"><li>Printer on</li><li>Please contact METTLER TOLEDO service</li></ul>
Blue Bluetooth display flashes continuously	<ul style="list-style-type: none"><li>No connection</li><li>Bluetooth module faulty</li></ul>	<ul style="list-style-type: none"><li>Establish connection</li><li>Reduce distance</li><li>Please contact METTLER TOLEDO service</li></ul>

## **7 Consumables and Spare Parts**

	<b>Type / Order number</b>	<b>Standard</b>
Paper roll, set of 5	00072456	2 piece
Paper roll, adhesive, set of 3	11600388	1 piece
Ribbon cartridge, black, minimum order quantity: 2	00065975	1 piece
Spare printing unit (may be changed only by METTLER TOLEDO service!)	00600133	–
Power line fuse 230 V (T100L250V) or 115 V (T200L250V) (may be changed only by METTLER TOLEDO service!)		commercially available

## 8 Technical Data

### Printer

Printing unit	Dot matrix printer 5 × 9 dots, 24-character line length, Epson IBM character table No. 4
Printing speed	1 line per second
Ribbon cartridge	Exchangeable, black
Paper roll	Normal paper, 58 × dia. 51 mm, integrated in housing, commercial size

### Printer functions

Printing, lightfast and thermally stable (GLP, GMP, ISO 9001), line feed, self-test

### Data transmission

Wireless Bluetooth® technology  
Max. range of 10 m

### Designed test conditions

Power line voltage, frequency	115 V or 230 V, admissible fluctuations –20% +15%, 50/60 Hz
Power consumption	140 mA or 70 mA
Range of application	For use only in closed interior rooms
Altitude	4000 m
Temperature range	+5 °C to +40 °C
Humidity	Max. relative humidity 80% for temperatures to max. 31 °C, linear decreasing to 50% at 40 °C
Installation category	II
Pollution degree	2
Dimensions / Weight	W × D × H = 157 × 210 × 85 mm / net 1.5 kg (incl. paper roll)

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© Mettler-Toledo AG 11/2013
### **Mettler-Toledo AG Laboratory & Weighing Technologies**