
Operating Instructions
Bedienungsanleitung
Mode d'emploi
Instrucciones de manejo

METTLER TOLEDO

PE200
PE400
PE2000
PE4000

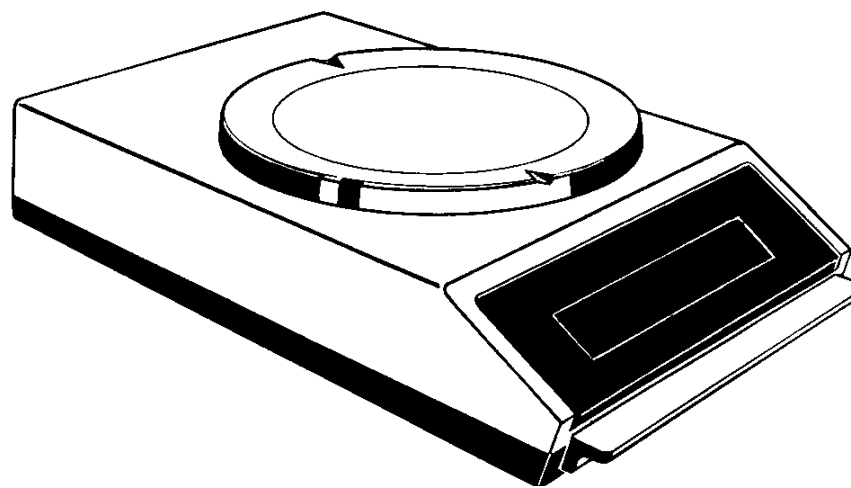OPERATING INSTRUCTIONS

Page 2

English

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Seite 20

Deutsch

MODE D'EMPLOI

Page 38

Français

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
Página 56

Español

AVAILABLE MODELS

These Operating Instructions apply to all models of the METTLER TOLEDO PE Series introduced on these pages. The operation of the balances is virtually identical. The differences are only apparent in their weighing range, readability and size of the weighing pan.

PE balances are equipped with METTLER TOLEDO DeltaDisplay, a special display mode for weighing in.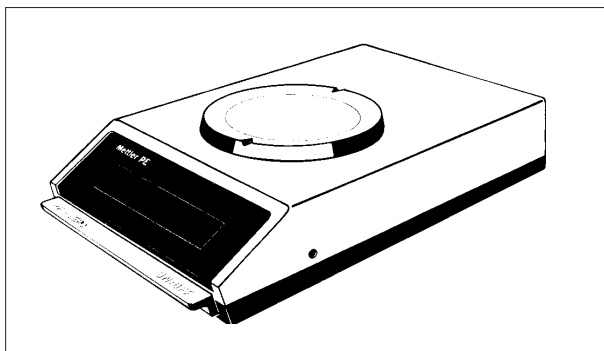


PE200

Weighing range 210 g
Readability 0.01 g

PE400

Weighing range 410 g
Readability 0.01 g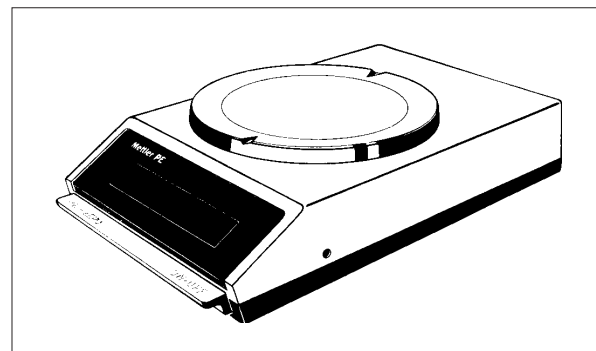


PE2000

Weighing range 2100 g
Readability 0.1 g

PE4000

Weighing range 4100 g
Readability 0.1 g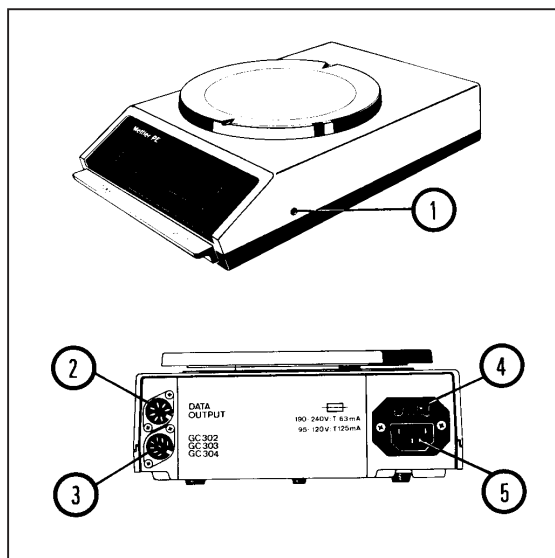

Operating components and connections

- 1 Calibration screw
- 2 Connection socket for Data Output (Option 05)
- 3 Connection socket for GC302, GC303 and GC304
Application Input Units
- 4 Fuse holder with back-up fuse
- 5 Power cable receptacle

TABLE OF CONTENTS

Page

Preparation

- How to check the voltage setting 4
- How to position the weighing pan and set up the balance 5

Operation

- How to switch the balance on or off 6
- How to calibrate the balance 7
- How to make a weight determination or tare the balance 8
- How to weigh in with DeltaDisplay 9
- How to weigh out and make weight comparisons 10

Additional features

- Unstable or restless weighing objects; weighing below the balance 11
- Application Input Units; 05 Data Output 12

Peripheral instruments (for 05 Data Output)

13

Accessories

14

Care and Maintenance

- How to replace the in-use cover and the fuse 15
- Maintenance and calibration 16

What if ...

17


Technical specifications

18

PREPARATION: How to check and correct the preselected voltage

The voltage selected for the balance must agree with the power line voltage. Before placing the balance in operating for the first time, check and if necessary, correct the voltage setting.


The factory setting is indicated on a label at power cable receptacle (5). If the indicated voltage agrees with that of the local power supply line, this page can be disregarded. If the voltage does not agree with the local power supply voltage, or if the label is missing, open the balance housing, check the position of the voltage selector and change it, if necessary. To do this, proceed as follows: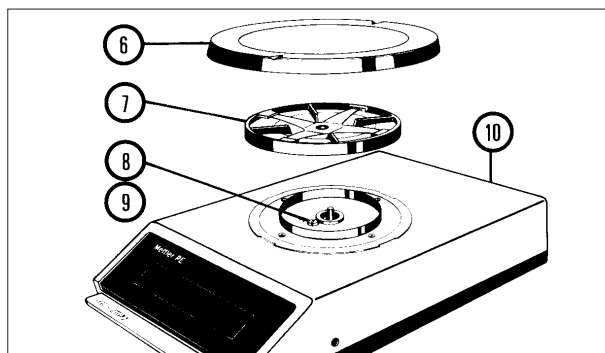

Opening the balance housing

Make sure the power cord is not connected

- Remove weighing pan (6) and pan support (7), if they are already installed.
- Remove screw (8) and put aside toothed ring and washer (9).
- Carefully lift off upper part of housing (10) together with in-use cover, if the latter is installed.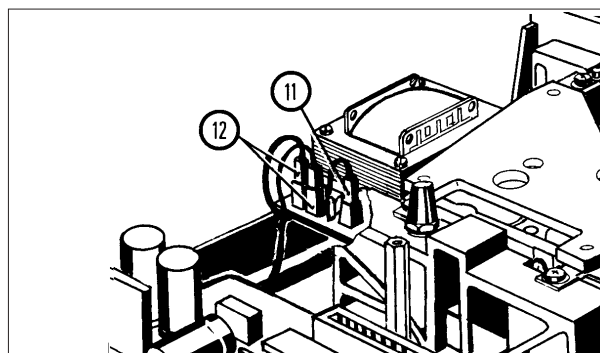


Voltage selector next to weighing cell

- Check to make sure that voltage selector plug (11) is attached to the pin which designated your local power supply voltage.
- If necessary, disconnect voltage selector plug (11) and attach it to the pin designated with your local power supply voltage.
- Admissible operating voltages:
 - 95 V, 105 V { insert
 - 110 V, 120 V { 125 mA microfuse
 - 190 V, 210 V { insert
 - 220 V, 240 V { 63 mA microfuse

Note: when changing the power supply voltage setting from 95...120 V to 190...240 V (or vice versa), make sure that the microfuse is also changed.

- The two connectors (12) must not be interchanged.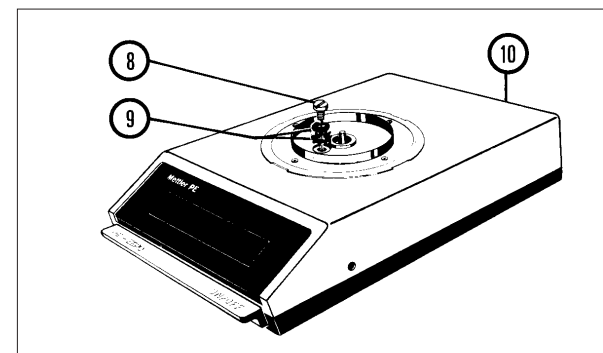




Closing the balance housing

- Carefully place upper part of housing (10) from top onto balance.
- First position toothed ring, then regular washer (9).
- Screw in and tighten screw (8).

PREPARATION: How to install the weighing pan and set up the balance

Installing the pan support and weighing pan is always done in the same manner, regardless of whether the in-use cover is installed or not.

Every PE balance provides reliable results, even under less than favorable environmental conditions. However, it is best to select a location that is stable, and not exposed to sunshine or to drafts (fast, stable display).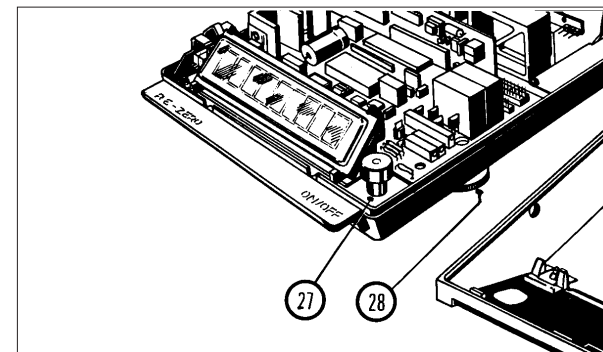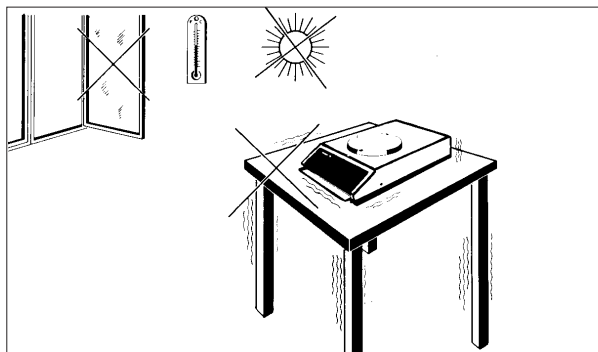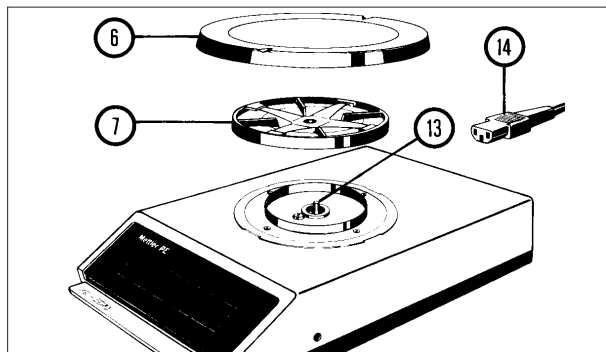

Installing the weighing pan

- Place pan support (7) on conical peg (13).
- Place weighing pan (6) on pan support.
- Connect power cable (14).

Location

- Solid, vibration-free support.
- No excessive temperature fluctuations.
- Avoid exposure to direct sunlight.
- Avoid drafts.

If it is absolutely necessary to work under less favorable conditions, consult section "ADDITIONAL FEATURES: Switching the integration time".


Level indicator field installation kit

This level indicator field installation kit was developed for PE balances which are moved frequently to new locations. With the level indicator, the PE balance can be leveled at every new location. Its efficiency is therefore increased. The field installation kit contains 2 foot screws (28) and 1 level indicator (27). METTLER TOLEDO Service can also be called on for this installation. Order No. see page 14.

OPERATION: How to switch the balance on or off

If you use your PE balance several times a day, it is best to leave it switched on all day long. This way, a constant operating temperature can be reached and weighing accuracy is increased.

It is also recommended that the balance be switched on some time before you begin to work with it (warm-up time).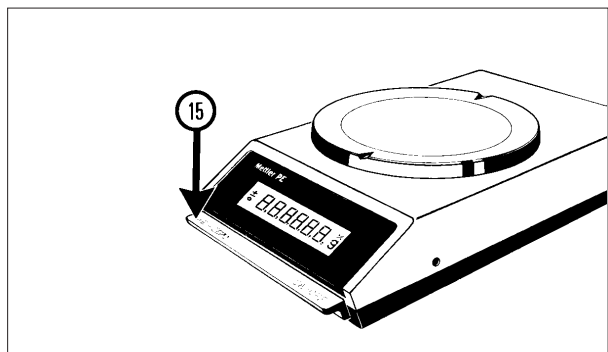


Switching the balance on

- Briefly press single control bar (15): all display elements light up for a few seconds.

± 888888.8 g

This will allow you to make a function check of the display. Then the zero display lights up with as many decimal places as are provided by the accuracy of your balance model,


- i.e. 0.00 g for the PE200, PE400
- 0.0 g for the PE2000, PE4000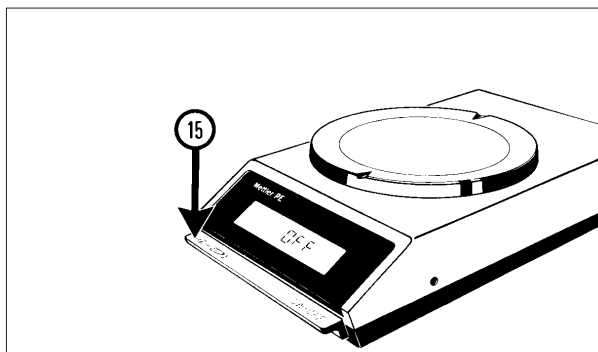

Display indicates OFF

If the power supply is interrupted during operation, the display will indicate "OFF" as soon as power is restored.

At that time, you must briefly press single control bar (15).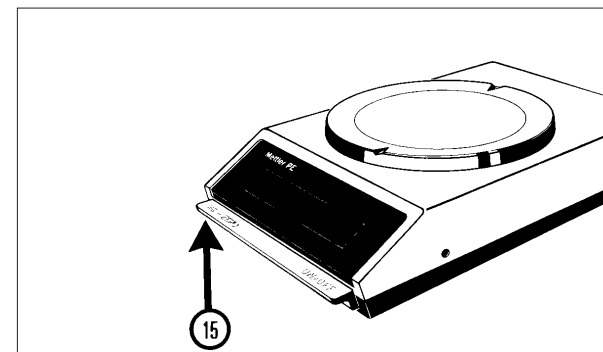


Switching the balance off

- Lift single control bar (15).

OPERATION: How to calibrate the balance

Your balance must be calibrated (i.e., set to the correct weight indication) every time its location is changed; this is the only way to ensure accurate weighings. Before being calibrated, the balance must be left switched on for at least 30 minutes (warm-up time).

Depending on your balance model, you will need certain test weights for calibration. These are listed under "ACCESSORIES, Optional equipment" for each balance model. The test weights must always be handled carefully; they should not be picked up with your bare hands.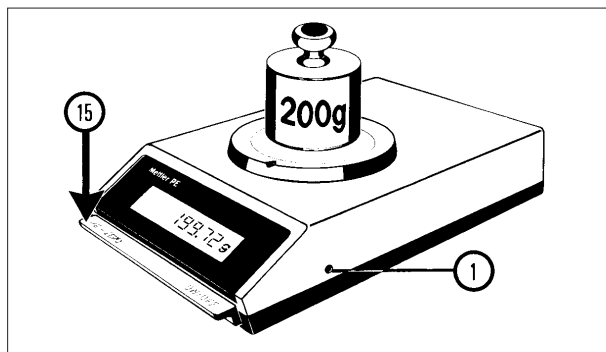


Checking the calibration

- Press single control bar (15): display will indicate zero.
- Place test weight on pan, read indication on display.

If the result displayed by the balance corresponds to the weight indicated on the test weight down to the last decimal, your balance is O.K. If not, it must be calibrated.

Using calibration screw (1), adjust calibration as described in the adjacent column.

Recheck calibration after every correction (see column at right).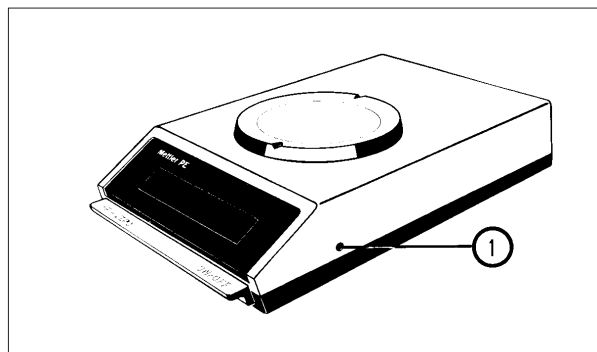


Adjusting the calibration

- Place test weight on pan, press single control bar.
- Remove test weight from pan.
- Turn calibration screw (1):
 - Clockwise, if the indicated weight value is larger than the test weight;
 - Counterclockwise, if the indicated weight value is smaller than the test weight.

With each turn of the screw, the calibration is changed by about:

0.16 g for the PE200, PE400 (test weight 200 g),
1.6 g for the PE2000, PE4000 (test weight 2000 g).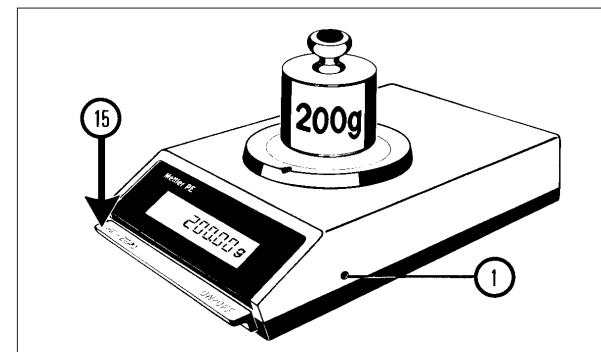

Rechecking the calibration

After every correction made with calibration screw (1), the calibration must be rechecked:

- Press single control bar (15), display will indicate zero.
- Place test weight on pan and read result.

If the result now displayed by the balance corresponds to the weight indicated on the test weight down to the last decimal, your balance is ready for use.

If not, continue adjusting calibration screw (1) until the balance is correctly calibrated.

OPERATION: How to make a weight determination or tare the balance


In PE balances, the readability is the same throughout the entire weighing range.

Automatic rounding of the last digit increases the accuracy of your PE balance even more.

Taring means that the display is reset to zero while a weight (such as a container) is on the weighing pan.

This way, the balance will automatically subtract the weight of the container from the total weight when an object or substance is weighed in. In that case, the balance only displays the weight of the weighed-in object.


However, the combined weight of container and the object may not exceed the overall weighing range.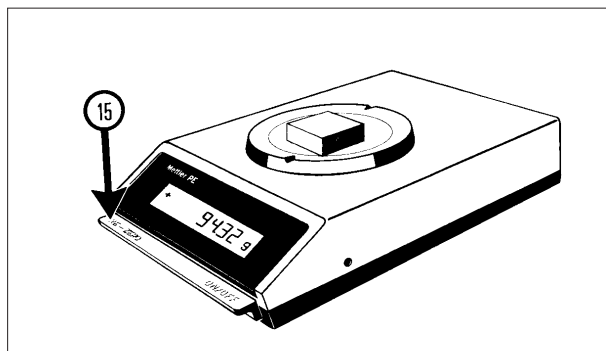

Weighing

- Briefly press down single control bar (15) to set the display to zero.
- Place weighing object on pan.
- Read weight display.

If the weighing range is exceeded, the display is blanked out except for the upper horizontal segments of the numbers. The balance thus indicates that it is overloaded.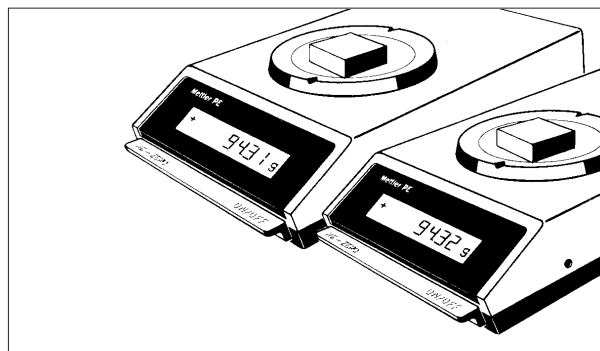


Display accuracy (result rounding)

Your balance always measures one more decimal place than it actually displays. This last, invisible digit is then rounded off according to the 4/5 principle.

Example:

The balance measures 94.314 g,
but displays 94.31 g

or
the balance measures 94.315 g,
but displays 94.32 g.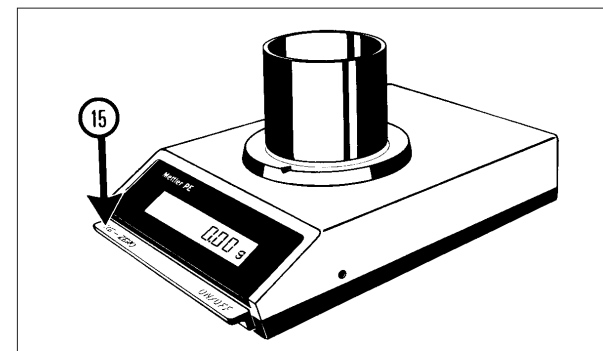

Taring

- Place container on weighing pan: balance will indicate its weight.
- Briefly press down single control bar (15), thus taring the balance: display now indicates zero.


The entire weighing range (minus the weight of the container) is now available for weighing in.

If DeltaDisplay is switched on (see next page), and if the balance is tared before achieving stability, the entire display is blanked out until stability is reached; only then will zero appear.

OPERATION: How to weigh in


Weighing-in means to fill loose or liquid objects into a container on the pan until a desired weight (target weight) is reached.

All PE balances are equipped with DeltaDisplay which considerably simplifies weighing in. To help you follow the weight changes, the display sequence speeds up automatically and the last decimal place of the display is blanked out. A stability detector is switched on along with the DeltaDisplay and it blocks the data output until the weighing result is stable enough. DeltaDisplay and stability detector can be switched off from within the balance (see also "ADDITIONAL FEATURES OF PE BALANCES").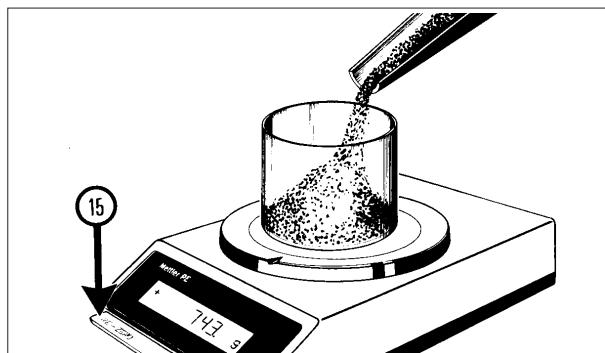

Weighing-in

- Place container on pan.
- Tare (15): Balance will now indicate zero.
- Fill in weighing objects until desired target weight is reached.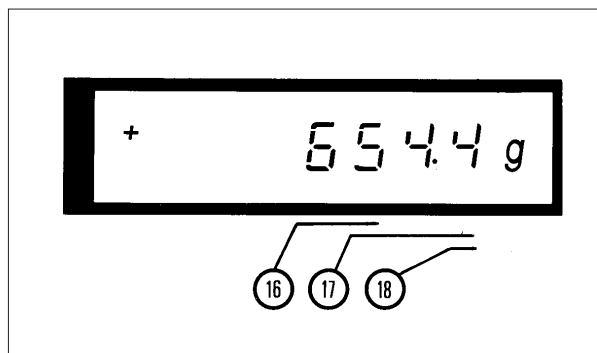


Reading the result

While filling in at a steady and fairly rapid pace until you reach the vicinity of the target weight, watch only the first two digits (16).

During fine weighing-in to the actual target weight, watch primarily the digits at right (17).

When the weight increases rapidly, there is also an automatic increase in the display speed and the last digit (18) is temporarily blanked out. It reappears again when fine weighing-in is resumed.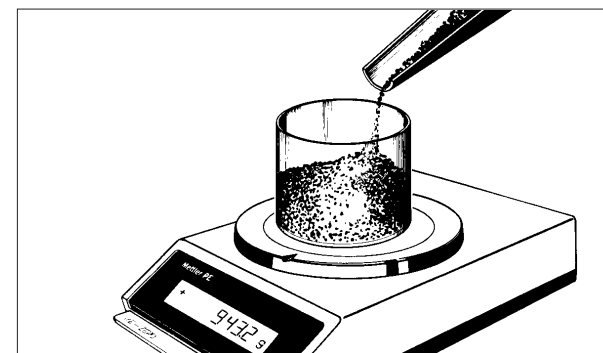

Weighing-in several components


If different objects are to be weighed in one after the other, the operator can tare after each component. This way, every new weighing can start from zero. This procedure can be continued until the combined weight of the container and of the weighing objects reaches the end of the total weighing range of the balance:

- PE200 210 g
- PE400 410 g
- PE2000 2100 g
- PE4000 4100 g

OPERATION: How to weigh out; how to read deviations from a reference weight


It is also possible to tare out a container filled with weighing objects. Thus the weighing-in process can be reversed and it is just as feasible to weigh out objects from a previously weighed full container.

Likewise, it is possible to tare a target weight (reference weight). This way, deviations from the reference weight can be read directly.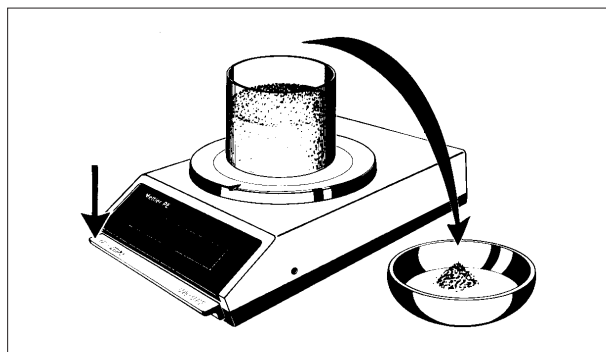

Weighing out


- Place container with weighing object on pan.
- Tare: Balance will now indicate zero.
- Remove material from container:
The weight of the removed material will now be displayed with the minus sign in front.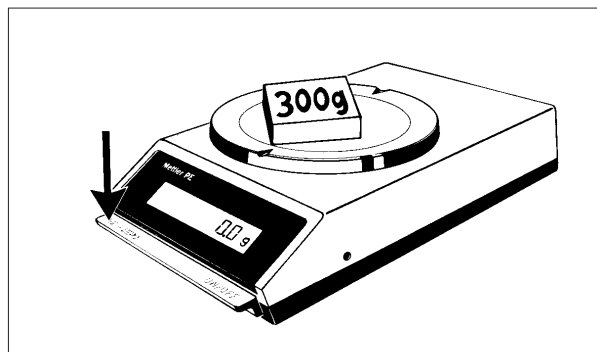

Entering a reference weight

- Place reference weight on pan (weight piece or sample).
- Tare: Balance will now indicate zero.
- Remove reference weight.
The reference weight will now appear in the display with the minus sign in front.
Example: -300.0 g

Deviations from this reference weight can now be read directly: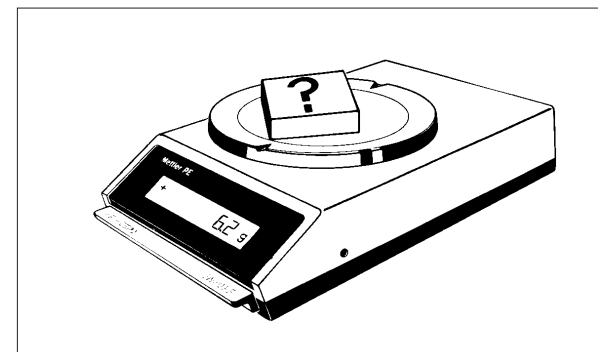

Determining a deviation

- Place object that is to be compared with reference weight on pan.


If the object is heavier, the actual deviation is indicated on the display with the positive sign in front, because:
 $-300.0 \text{ g} + 306.2 \text{ g} = +6.2 \text{ g}$

If the object is lighter, the actual deviation is indicated with the negative sign in front, because:
 $-300.0 \text{ g} + 294.2 \text{ g} = -5.8 \text{ g}$

ADDITIONAL FEATURES OF PE BALANCES: Restless weighing objects; weighing below the balance

All PE balances are shipped with the DeltaDisplay switched on. In case of rapid weight changes, the display changes are automatically increased and the last digit is temporarily blanked out. The DeltaDisplay can be switched off by an internal switch. In addition, the integration time can be doubled. In this case, display changes are only half as fast. This makes it possible to adapt PE balances to difficult conditions and to stabilize restless displays to the correct value.

All PE balances are equipped for weighing below the balance.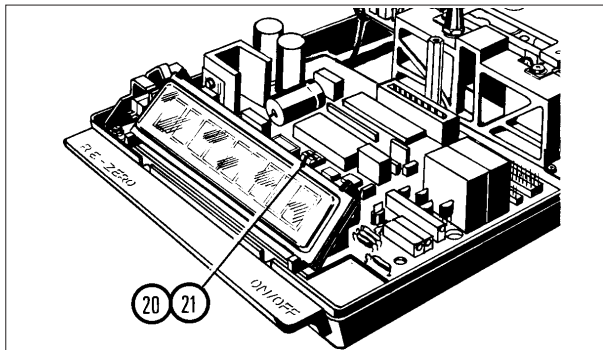


Switching DeltaDisplay and integration time

If restless objects have to be weighed (e.g., live animals) or if weighing must be made under conditions of external instability (i.e., drafts), the result can be stabilized by switching of DeltaDisplay and doubling the integration time.

This procedure can also be handled by METTLER TOLEDO Service.

- Disconnect power cable.
- Open balance housing (see page 4, "PREPARATION; How to check the voltage setting").

Switches (20) (21) for DeltaDisplay and integration time are located directly behind the display.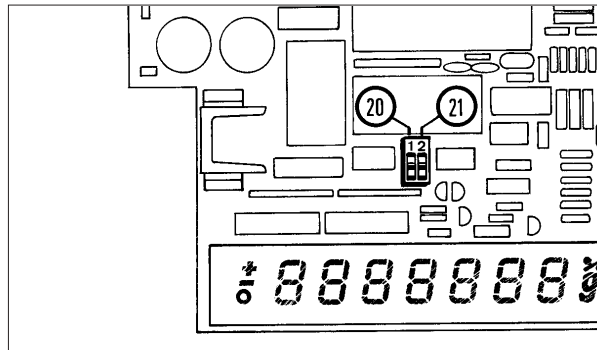


Slide switch (20) at left controls DeltaDisplay and stability detector.

- DeltaDisplay and SD on: Switch is at the stop nearer to the display.
- DeltaDisplay and SD off: Switch is at rear stop.

Slide switch (21) at right is used to adjust the integration time:

- Short integration time: Switch is at the stop nearer to the display.
- Long integration time: Switch is at rear stop.
- Close balance housing (see under "PREPARATION: How to check the voltage setting").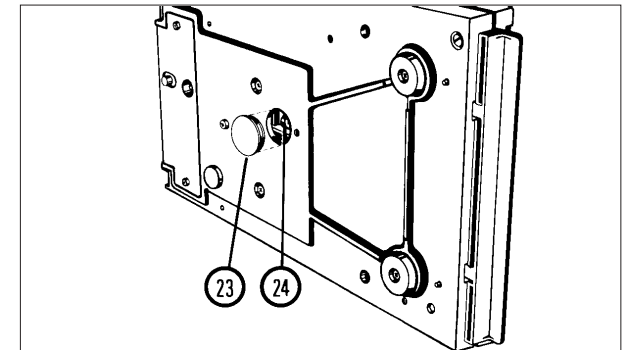

Weighing below the balance


Your PE balance is specially equipped to handle weight determinations that can not be made on the weighing pan. This is done by suspending the object from the weighing cell. To do this, the balance table must have an appropriate opening.

- Remove weighing pan and pan support, and tilt balance to its side.
- Remove cover (23).
- Using a holder, suspend weighing object from hook (24).
- Set balance right side up and install weighing pan and pan support.
- With holder attached, tare balance.

ADDITIONAL FEATURES OF PE BALANCES: Application Input Units; 05 Data Output

Some very useful accessories for the PE200/400/2000/4000 are the GC302, GC303 and GC304 Application Input Units. They make it possible to quickly and conveniently perform a variety of special weighing procedures and functions.

A data output (option 05) can be delivered along with every PE200/400/2000/4000 and easily installed by the user. The necessary wires and connectors are already installed at the factory. In the next two pages we show you some of the METTLER TOLEDO peripheral instruments and accessories which make it possible to use PE balances for special applications.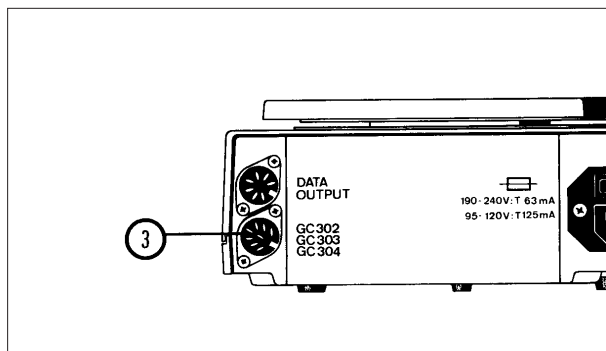

Connection for Application Input Units


Every PE200/400/2000/4000 is equipped with a connection socket (3) for GC Application Input Units.

Application Input Units

These units are suitable for the following applications:

GC302 Application Input Unit

- Piece counting
- Percent weighing
- Net total weighing
- Animal weighing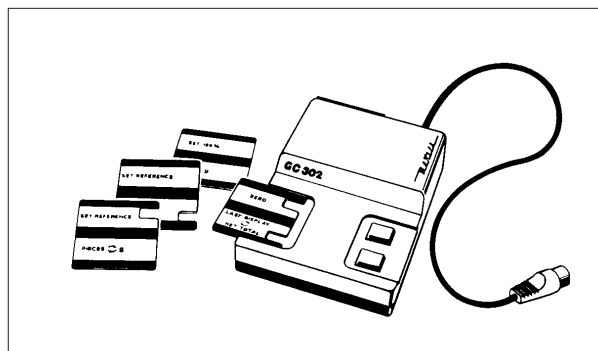


GC303 Application Input Unit

- Piece counting with weight display in:
pounds/grams (lb/g)
ounces/grams (oz/g)

GC304 Application Input Unit

- Weighing in grams and in the following non-metric units:
 - Pounds (lb)
 - Ounces (oz)
 - Troy-ounces (ozt)
 - Pennyweights (dwt)
 - Grains (GN)
 - Tael (tl)
 - Carats (C.M.)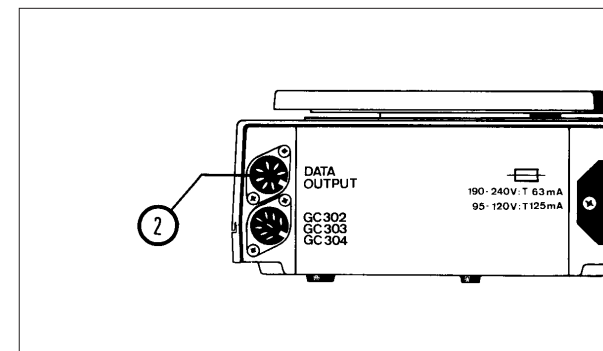

05 Data Output

The following peripheral instruments can be connected to the 05 Data Output (2):

- GA40 Thermal Printer
- GC45 Remote Display
- GC47 Digital/Analog Converter


(These instruments are discussed in greater detail on the following page).

You can also retrofit your balance by using an "Option 05 Field Installation Kit" (Order No. 47660).

PERIPHERAL INSTRUMENTS: Data output and data preparation

The three peripheral instruments introduced on this page can be connected to any PE200/400/2000/4000 balance equipped with a 05 Data Output (see preceding page).

Several peripheral instruments can be connected at the same time.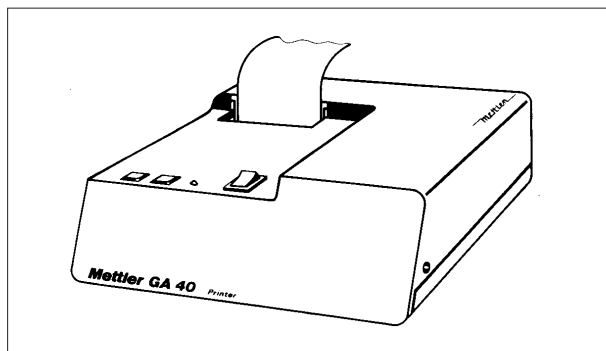


GA40 Thermal Printer

For the printed recording of weighing data.
Maintenance-free thermal printer (dot printer, no printing ribbon), compact, fast and low noise.

Nonfading paper tape, not sensitive to mechanical imprinting, rolls are 30 meter in length each and will accommodate about 5000 lines of print. Width of paper tape 57 mm (2 1/4 inch).

Separate power supply:
110/130/220/240 V

Dimension (W x D x H):
185 x 200 x 60 mm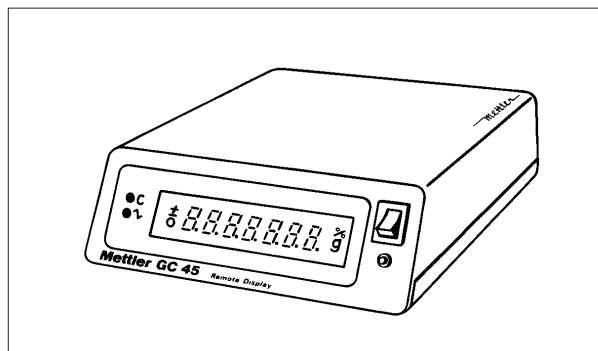


GC45 Remote Display

For the visual monitoring of weighing results away from the weighing station.

Display is completely identical with that of the balance itself.

Separate power supply:
110/130/220/240 V

Dimension (W x D x H):
185 x 200 x 60 mm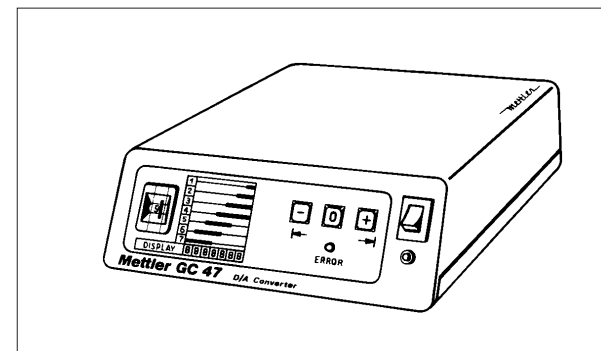

GC47 Digital/Analog Converter

For the conditioning of (digital) weighing data generated by the balance for entry into analog instruments such as line printers.

Two outputs: ± 10 V; ± 1 V


Multiswitch for decade selection (in group of three).

Separate power supply:
110/130/220/240 V

Dimension (W x D x H):
185 x 200 x 60 mm

ACCESSORIES


Optional equipment: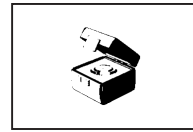

Calibration weights
certified according to OIML
200 g for PE200/400
2000 g for PE2000/4000


Order No.

42053
42065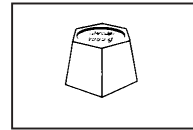


Calibration weights, non-certified
200 g for PE200, PE400
2000 g for PE2000, PE4000

47698
47699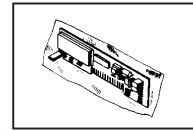


Field Installation Kit
for 05 Data Output

47660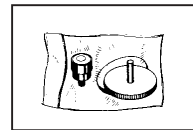


Field Installation Kit
for level indicator

47659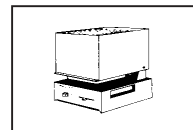


Infrared Dryer
110/220 V, 140 VA

LP15B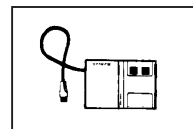


Application Input Unit

GC302
GC303
GC304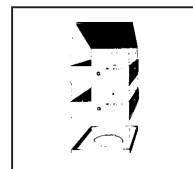

Lid 46433
Element 46432
Base 43256 } Plastic draft
shield with two
elements
(can only be used with PE200/400)


43260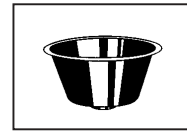

Animal weighing pan
(for PE2000/4000 only)

Order No.


47644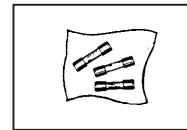

Microfuses (set of 3)
125 mA, slow-acting, for 95/120 V
63 mA, slow-acting, for 190/240 V


26172
46328

Standard equipment: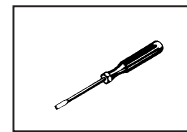


Screwdriver

50279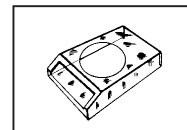


In-use cover

47465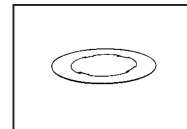

Retaining ring for in-use cover

41156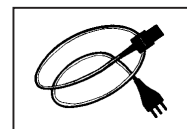

Power cable

Neutral
Switzerland
Germany
USA

87576
87920
87925
88668


CARE AND MAINTENANCE: How to replace the in-use cover and the fuse

To protect the balance housing against soiling, damage and corrosive chemicals, the balance is delivered with the in-use cover already in place. This cover is made of stiff plastic so that it cannot be folded.

The in-use cover does in no way hinder your working with the balance.


A back-up fuse is built into the fuse holder, in case your microfuse becomes defective.

If the microfuse has to be replaced repeatedly without apparent reason, call METTLER TOLEDO Service.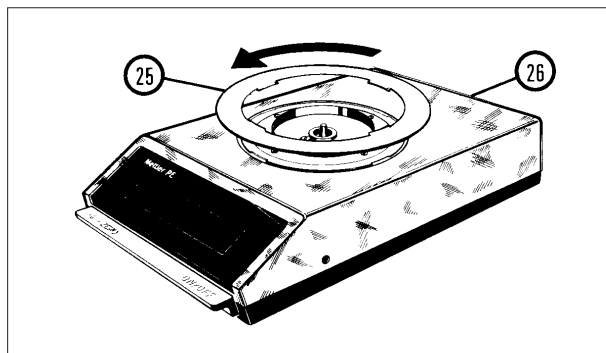


Removing the in-use cover

- Remove weighing pan and pan support.
- Turn retaining ring (25) counterclockwise until it is released and can be lifted off.
- Remove in-use cover (26).
- Install pan support and weighing pan.
(The retaining ring should only be installed when the in-use cover is in place).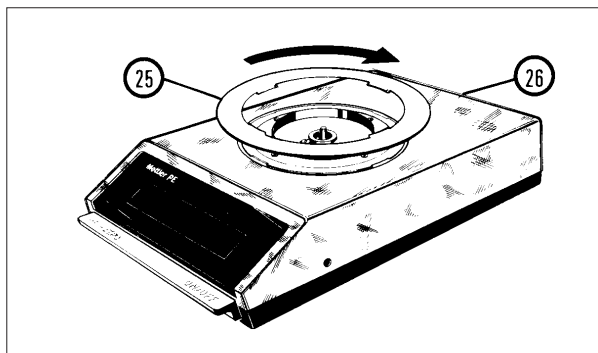

Installing the in-use cover

- Remove weighing pan and pan support.
- Place in-use cover (26) on balance.
- Position and engage retaining ring (25), then turn clockwise to stop position.
- Install pan support and weighing pan.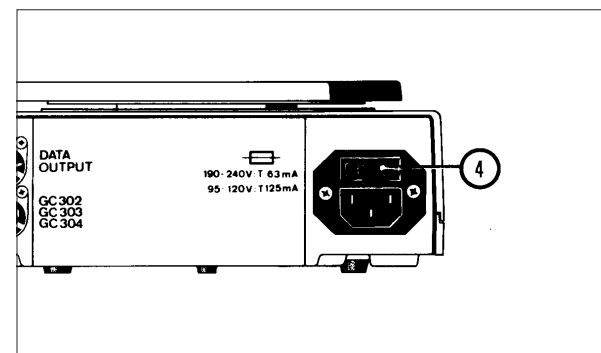


Replacing the microfuse

- Disconnect power cable.
- Using a screwdriver, remove fuse holder (4).
- Remove defective fuse and insert new fuse (back-up fuse from fuse holder).
Fuse rating:
125 mA, slow-acting, for 95/120 V
63 mA, slow-acting, for 190/240 V
- Insert fuse holder.
- Plug in power cable.

CARE AND MAINTENANCE: Upkeep and calibration

Your METTLER TOLEDO PE balance requires very little maintenance, but it should be cleaned and calibrated at regular intervals.


Before checking the calibration, the balance must have been switched on for at least 30 minutes (warm-up time).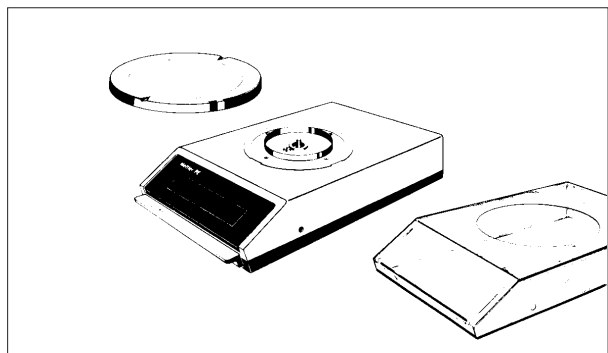

Cleaning

Weighing pan, in-use cover and, if necessary, the balance housing should be cleaned whenever the need arises.

As much as possible, avoid using strong solvents.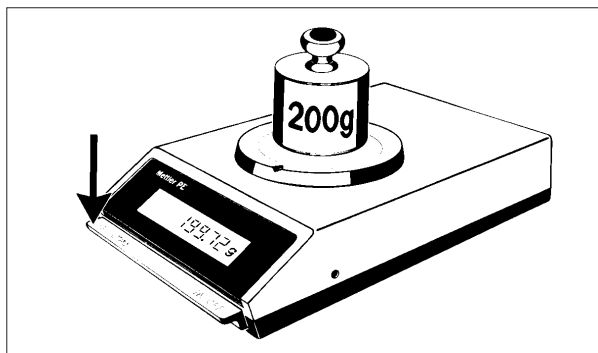


Calibration

If your balance is to deliver accurate weighing results, it must be properly calibrated.

Check the calibration of your balance at regular intervals; in case of constant operation, at least once a month.

The test weight to be used with your balance model is listed under "ACCESSORIES, Optional equipment".

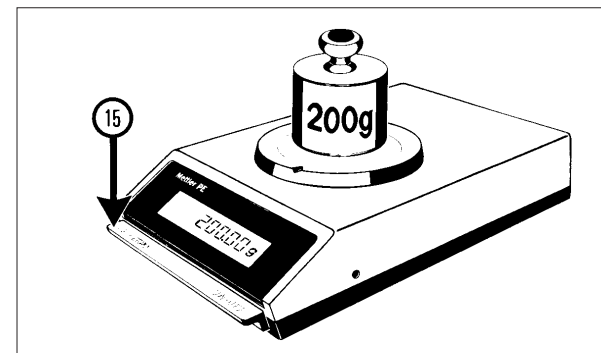Rechecking the calibration

- Press single control bar (15):
The display indicates zero.
- Place test weight on pan and read the result.

If the result now displayed by the balance corresponds to the weight indicated on the test weight down to the last decimal, your balance is ready for use. If not, it has to be calibrated.

For instructions on how to do it, see "OPERATION: How to calibrate the balance" (page 7).

WHAT IF ...

... the entire display is blanked out?

... only the upper segments of the display light up?

... only the lower segments of the display light up?

... OFF is displayed?

... the weighing result is unstable?

... the weighing result is obviously wrong?

Then ...

- Balance is not switched on.
- Power cable is not plugged in.
- No power.
- Microfuse is defective (see "CARE AND MAINTENANCE: Replacing the microfuse").
If this occurs again: Check voltage selector setting and fuse rating.
If both are correct, call METTLER TOLEDO Service.
- Weight on pan is too heavy (weighing range is exceeded).
- Balance is defective (call METTLER TOLEDO Service).
- Pan support is not installed.
- Weighing pan is not installed.
- In-use cover is touching pan support (retaining ring of in-use cover is not installed).
- Weighing range was exceeded before balance was unloaded (overload was not obvious because it is composed of tare and weighing object).
Correct defect by taring after the balance is unloaded.
- Power was out: Check the two power cable connectors for tight fit.
Again switch on balance by pressing control bar.
- Control bar was not pressed down all the way when power was switched on.
- Draft at the weighing station.
(PE200, PE400: Use draft shield. See "ACCESSORIES: Optional equipment").
- Balance table is not stable.
- Weighing object is restless, e.g., live animals (see "ADDITIONAL FEATURES").
- Voltage selector is set to an incorrect voltage (see "PREPARATION: How to correct the selected voltage setting").
- Balance is not leveled.
- Balance is not properly calibrated (see "OPERATION: How to calibrate the balance").
- In-use cover touches the pan support (install retaining ring).
- Balance was not tared before weighing was made (single control bar had not been pressed).
- Weighing object might be touching the draft shield.

TECHNICAL SPECIFICATIONS

	<u>PE200</u>	<u>PE400</u>	<u>PE2000</u>	<u>PE4000</u>
Weighing range	210 g	410 g	2100 g	4100 g
Readability	0.01 g	0.01 g	0.1 g	0.1 g
Tare range (by subtraction)		210 g	410 g	2100 g 4100 g
Admissible ambient temperature (during operation)	10...+ 40 °C	10...+ 40 °C	10...+ 40 °C	10...+ 40 °C
Sensitivity drift (10...30 °C)	$\pm 2 \cdot 10^{-5} / ^\circ\text{C}$	$\pm 1.2 \cdot 10^{-5} / ^\circ\text{C}$	$\pm 2 \cdot 10^{-5} / ^\circ\text{C}$	$\pm 1.2 \cdot 10^{-5} / ^\circ\text{C}$
Result deviation (with balance inclined 1 : 1000)	± 0.02 g	± 0.02 g	± 0.2 g	± 0.2 g
Reproducibility	0.005 g	0.005 g	0.05 g	0.05 g
Linearity	± 0.015 g	± 0.015 g	± 0.15 g	± 0.15 g
Stabilization time (typical)		2 s		
Integration time		1.25 s / 2.5 s*		
Display sequence: DeltaDisplay off		0.6 s / 1.2 s*		* When integration time is doubled
DeltaDisplay on		0.15 s / 0.3 s*		
Stability control (sensitivity invariable)		ON / OFF		
Power supply:				
– Adjustable voltage settings		95V, 105V / 110V, 120V / 190V, 210V / 220V, 240V		
– Admissible voltage fluctuations		+ 10 %		
– Frequency		50...60 Hz		
– Power consumption		about 10 VA		
Dimensions of housing (W x D x H)		194 x 310 x 65 mm		
Weight	3.7 kg		3.8 kg	
Weighing pan (chrom-nickel steel)	dia. 130 mm		dia. 170 mm	

METTLER TOLEDO 05 Data Output Combined current loop and voltage interface.

Galvanically separated, unidirectional 2-wire interface. Characters: alphanumeric. Character code: ASCII/ISO, 7-bit, even parity.

Transfer speed: 2400 baud. Connector socket: 7-pole, type Hirschmann Mab 7S.

03 Mode (for METTLER TOLEDO peripheral instruments) Loop current: 20 mA (passive)

CL Mode (for METTLER TOLEDO CL instruments as well as for non-METTLER TOLEDO instrument with active loop current input)

RS232 Mode (for non-METTLER TOLEDO instruments) Voltage interface according to RS232C/V24

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© Mettler-Toledo GmbH 2001 701031 Printed in Switzerland 0111/2.33
Mettler-Toledo GmbH, Laboratory & Weighing Technologies, CH-8606 Greifensee, Switzerland