
Operating instructions
Bedienungsanleitung
Mode d'emploi
Instrucciones de manejo
Istruzioni d'uso

METTLER TOLEDO

Printer METTLER TOLEDO Sprinter 1
Drucker METTLER TOLEDO Sprinter 1
Imprimante METTLER TOLEDO Sprinter 1
Impresora METTLER TOLEDO Sprinter 1
Stampante METTLER TOLEDO Sprinter 1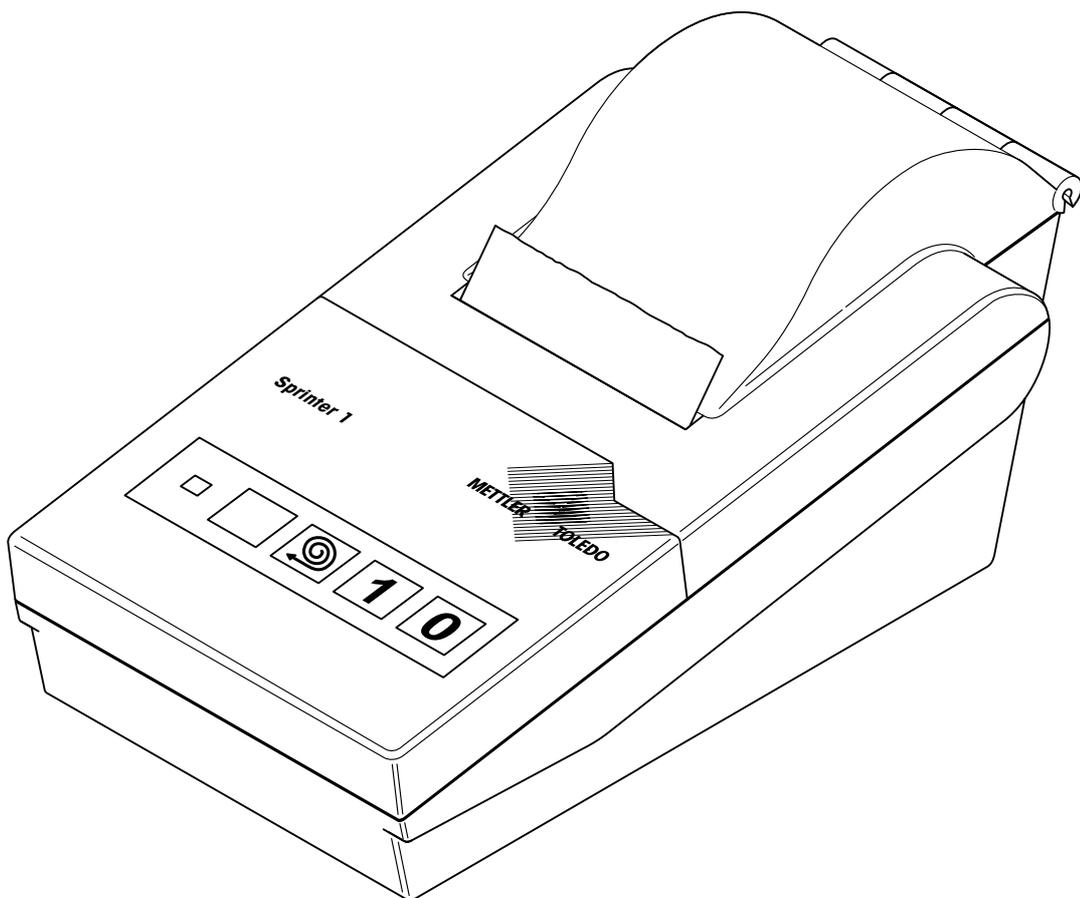

1. Safety

- Read through the operating instructions carefully.
- Check the standard equipment: Printer with ribbon and roll of paper, plug-in power supply unit, operating instructions.
- The printer must not be operated in wet areas or in a hazardous environment.
- Use only the plug-in power supply unit supplied and ensure that the voltage value printed on it matches your local line voltage. If this is not the case, on no account put the printer into operation.
- Use only optional equipment and expendables supplied by METTLER TOLEDO.
- As the printer is a precision unit, take appropriate care when handling it.
- Never open the printer, it contains no parts which can be repaired or changed by the user.

2. Keys and LED

Keys 1 and 0: On/off switching.

Feed key (key with paper roll symbol): Advance of the printer paper.

Programming key (key without a symbol): Configuration of the printer (see section 5).

Green LED: Printer ready for operation.

3. Preparation of the printer

- Insert paper roll in the printer. Push leading edge of paper as far as it will go into the slot below the ribbon cassette. Switch printer on with the 1 key and press feed key. Lead paper through slot in the cover and close cover.
- Connect power supply unit to the printer and the power supply. When putting into operation for the first time, the built-in battery must be charged via the power supply unit (approx. 14 hours).
- Connect interface cable to the printer and the balance.
- Configure the interface of the balance as follows (see balance instructions):
 - Operating mode: Printer *
 - Protocol: XON/XOFF *
 - Data bits/parity: 7/even *
 - Data transmission rate: 2400 baud ** = factory setting

4. Printer test

- With the printer switched off, press and hold the feed key and switch the printer on with the 1 key. The entire character set and the preselected settings will be printed out.
- Switch printer off with the 0 key to end the test.

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5. Configuring printer

- With the printer switched off, press and hold the programming key and switch the printer on with the 1 key.
- Use the feed key to select the parameter to be changed. After every keystroke, the current setting of the corresponding parameter will be printed out.
- Press the programming key to change the current setting of the selected parameter. Each time the key is pressed, the setting is changed and printed out. **Recommendation:** Work with the factory settings (see printout opposite).
- Use the feed key to select the next parameter or by pressing the programming key and the feed key simultaneously store the settings and exit the configuration mode (all settings will be printed out).

6. Changing the ribbon

- Open printer cover and press the ribbon cassette at the point marked "PUSH".
- Insert new cassette and press downward. Turn knurled wheel on cassette in the direction of the arrow to tension the ribbon. Close printer cover (pull paper through slot in the cover).

7. Cleaning

- Wipe the exterior of the printer with a moist cloth and a mild commercial household cleaning agent and dry it immediately.

8. Spare parts, optional equipment, expendables

Designation	Number
Power supply Euro (230 V)	21253400
Power supply UK (240 V)	21253401
Data cable	21253677
Paper roll (set of 5)	72456
Ribbon	65975

9. Data of the power supply

Power supply Euro:	
Primary:	230 V AC/50 Hz
Secondary:	9 VDC/1.3 A
Power supply UK:	
Primary:	240 V AC/50 Hz
Secondary:	9 VDC/1.3 A

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Declaration of conformity

We hereby declare under our sole responsibility that the product **"Sprinter 1"** S/N: > **04B.2939.014** to which this declaration relates is in conformity with the following directives and standards:

Directive	Standard
73/23/EWG Low voltage	EN60950 Safety requirements
89/336/EWG Electromagnetic compatibility	EN55022 Radio interferences EN50082-1 Immunity
90/384/EWG for nonautomatic weighing instruments	EN45501 for nonautomatic weighing instruments

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Roland Schmäder, General Manager

Dr. Andreas Rinn, Quality Manager

To protect your METTLER TOLEDO product's future:

METTLER TOLEDO service assures you of quality, measuring accuracy and preservation of value of the METTLER TOLEDO products for years to come. Please send for full details of our attractive terms of service. Thank you.

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