
**Operating Instructions**  
**Bedienungsanleitung**  
**Mode d'emploi**  
**Istruzioni d'uso**  
**Instrucciones de manejo**

**METTLER TOLEDO**  
**NetEasy-M**

**METTLER TOLEDO**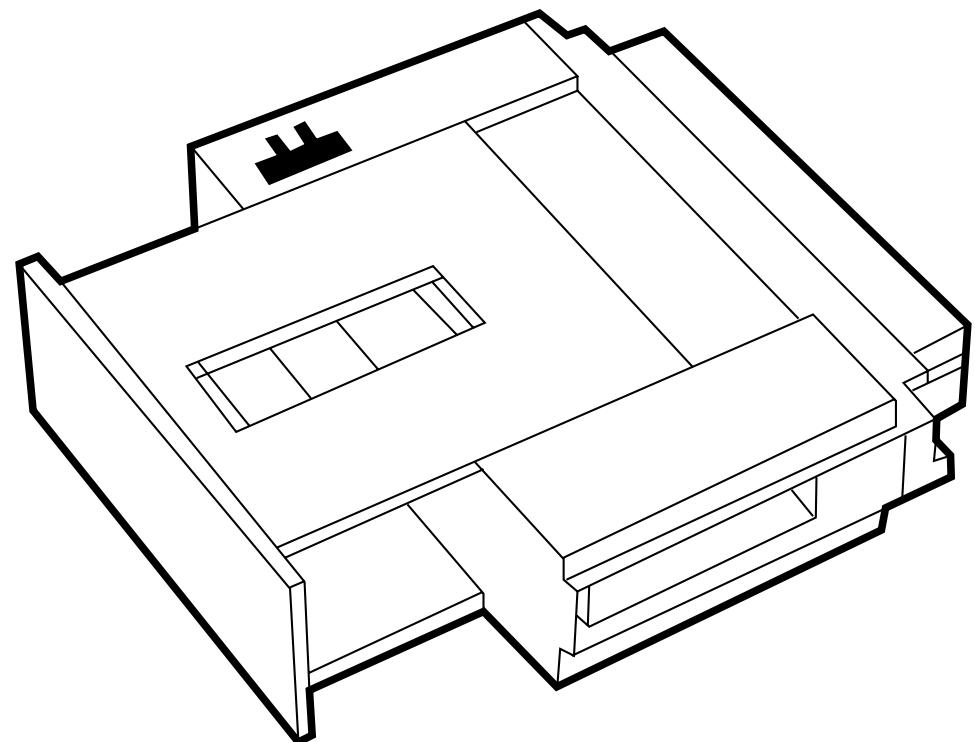

English Page 2

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
## General

NetEasy-M is a program cassette for SM scales, high-capacity PM scales or for AM/PM balances up to 6.1 kg with the GM303 Control Unit. It allows you to use your balance or scale for convenient net total weighing or totalization.

The **insertion of the program cassette** is described in the operating instructions of the balance or scale in the section “Configuration”. When NetEasy-M is inserted, the application “Net Total” is automatically activated after the balance or scale has been switched on.

## Configuration register with NetEasy-M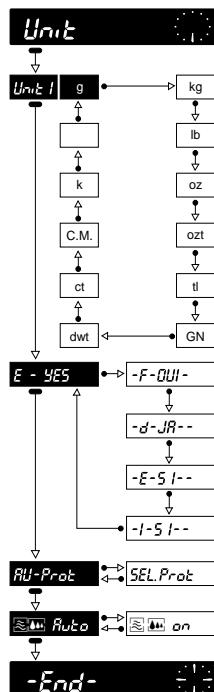

The configuration register with NetEasy-M is distinguished from the configuration register of the balance or scale with the standard program cassette only in the –Unit– sector. With NetEasy-M the following can be selected in the –Unit– sector.

Language	-E-YES-	English (default setting)
	-F-OUI-	French
	-d-JA-	German
	-E-SI-	Spanish
	-I-SI-	Italian


Printout	Au-Prot	automatic printout (default setting)
	SEL.Prot	printout only with the <b>PRINT</b> key

All further setting options are identical to those with the standard cassette, see operating instructions of the balance or scale.

### Notes

- If the configuration register can not be selected, check whether the jumper of the program cassette is in the position “unsecured” (see operating instructions of the balance or scale, section “Configuration”).
- When the balance or scale is reset to the default configuration (Reset), the language is not reset.

## Totalization

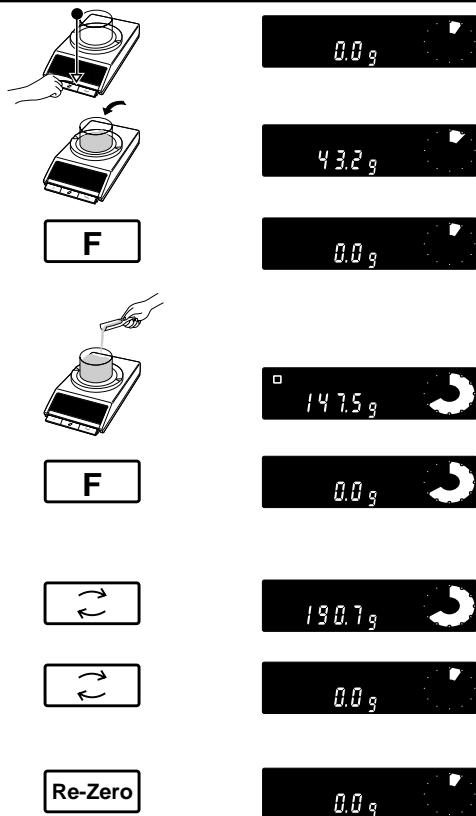NET TOTAL

-----  
Weight 43.2 g  
Weight 147.5 g  
Total 190.7 g  
-----

Load empty container, tare.

Add first component.

Store first component weight: press key **F**.  
This automatically resets the display to zero.

Add second (additional) component.

Store second (additional) component: press key **F**.  
This automatically resets the display to zero.

Switching between last display and current total (net total): press key **F**.

Recall of the last display: press key **F**.

Clear total (net total): press control bar (Re-Zero).

### Notes

- Each component weight is automatically printed out on storage if –Au-Prot– is configured with an attached printer.
- Printout of the net total weight: on display of the total (net total), press the **PRINT** key.
- If the total is displayed, the last component can be added up to the total target weight. The weight of the last component is not, however, listed when the net total is printed out.

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© Mettler-Toledo GmbH 2000 704713 Printed in Switzerland 0008/2.45
**Mettler-Toledo GmbH, Laboratory & Weighing Technologies, CH-8606 Greifensee, Switzerland**