CytoDirect

Automated cell counter

METTLER TOLEDO

1 Use & limits

The instrument is designed to be used in laboratories by trained staff and is intended to count the number of cells and assess their viability. The typical work environment for the instrument is a life-science laboratory, often a quality control laboratory (i.e., near-production) in academic and industrial settings. It can provide information on the cell count (CC) and cell viability (VIA) of various cell types in suspension. For more information, see the following link: www.mt.com/CytoDirect-Cell-List
The instrument is suitable for samples with the following characteristics:

- Compatible with the substances that the samples typically come in contact with.
- Free of air bubbles.

Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

Prohibitions on use

The instrument should not be used:

- For medical diagnostics.
- In potentially explosive environments.
- With hazardous substances (corrosive, flammable).
- With hot liquids (> 60 °C).

Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties. Mettler-Toledo GmbH assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. Mettler-Toledo GmbH assumes that the instrument owner provides the necessary protective gear.

2 Warnings

Safety notes

WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO AC/DC adapter designed for your instrument.
- 2 Keep all electrical cables and connections away from liquids and moisture.
- 3 Check the cables and the plugs for damage and replace damaged cables and plugs.

CAUTION

Slight injuries from exposure to laser radiation

A laser LED (Class I) is used as a light source in the instrument. If the light source is activated when no slide or slide adapter is inserted, looking directly into the laser light can cause slight injuries to the eyes.

- 1 Never look directly into the light source.
- 2 Discontinue the use of the instrument immediately.
- 3 Seek support for repair or replacement of the instrument.

NOTICE

Damage to the instrument or malfunction due to the use of unsuitable parts.

Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

Storage and Transport conditions

- Temperature: -10 °C to 50 °C.
- Relative humidity: 10% to 90% (non-condensing).

Instrument misuse

Make sure to adequately fill the sample carrier (at least up to the checkmark) before measuring to avoid the formation of air bubbles. Trapped air in the sample area will lead to incorrect measurement results. Clean the microscope sensors frequently to avoid error messages and artifacts in the images.

3 Notes & Instructions

Installation

Unpack the cell counter.

- Remove the cell counter from the protective packaging. Store the protective packaging for later transport over long distances.

- Check that you have received all parts listed in the scope of delivery. Inspect the parts visually for flaws or damage. If parts are missing or damaged, report it to your authorized METTLER TOLEDO organization.

Position the instrument.

- The instrument has been developed for indoor operation in a well-ventilated area.
- The following site requirements apply:
 - Ambient conditions within the limits specified in the technical data.
 - No powerful vibrations, direct sunlight, corrosive gas atmosphere, explosive atmosphere, powerful electric or magnetic fields.
- Procedure:
 - Place the instrument on a stable, flat, and fire-resistant surface.
 - Make sure that nothing blocks the air vent of the instrument.

First startup.

- Connect the charger and press the power button to start the instrument for the first time.

Charge the device.

- Recharging an empty battery takes approximately 3 hours.
- The cell counter can be switched on and used for measurements during the charging process.

WARNING

A document named „User Manual“ is available for this instrument. A QR code for the User Manual download is delivered with the instrument. The digital User Manual contains a description of the instrument and its use. Keep this document for future reference. Include this document if you transfer the instrument to other parties. Only use the instrument according to the User Manual. If you do not use the instrument according to this document or if the instrument is modified, the safety of the instrument may be impaired, and Mettler-Toledo GmbH assumes no liability.

The User Manual for CytoDirect™ can be downloaded by scanning the QR code found printed on the envelope contained in the device packaging or by visiting the following link: <http://www.mt.com/CytoDirectUserManual>

A hard copy of the User Manual of CytoDirect will be provided upon request. Please send your inquiries to your authorized METTLER TOLEDO organization. If electronic access to the User Manual cannot be ensured, the user is advised to print the document out before use.

Disposal and recycling

- This measuring device and its components contain electronic parts.
- EU Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) specifies that used devices should not be disposed of as general household waste, laboratory waste or hospital waste. Used systems must be collected separately, to increase the recycling rate and to reduce the impact on human health and on nature.
- The device is equipped with a LiPo battery. Batteries require special care in handling and disposal. Improper disposal may have adverse effects on human health and may harm the environment. This also applies to countries outside the EU, as per their specific requirements.
- At the end of its product life cycle, the device should be disposed of as electronic waste. Please contact your authorized METTLER TOLEDO organization. We will dispose of it in a professional and environmentally friendly manner. Disposal via municipal collection points is not permitted.
- The sample carriers used for measuring purposes must be disposed of after use in accordance with the customer's own laboratory regulations. If the sample carriers have been mixed with human or animal samples, special care is required when handling them. These samples may be contaminated with pathogenic germs or viruses and must be handled as potentially hazardous.

Links and Addresses

Hereby, Mettler Toledo declares that the radio equipment type in CytoDirect is in compliance with Directive 2014/53/EU. The full text of the EU declaration of conformity is available at the following internet address: <http://www.mt.com/CytoDirect-DoC>

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© 01/2024 METTLER TOLEDO. All rights reserved.