## Switching the instrument on/off

24.7°C ATC

(A)

M 0003

• A

Cal 1



## Changing the mode



Pressing MODE changes the mode: pH, mV, rel. mV.

## Performing a pH calibration

Switch to pH mode. Cal ኩ Change pH buffer group




Blinking A: mea-

Insert electrode




25.0°C MTC suring

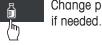

-2013 14:32 /A 25.0°C MTC ⑧/Ā 7.00 Value stable a 2.00, 4.01, 7.00, 9.21, 11.00 Call Exit End

Insert electrode into second pH buffer.

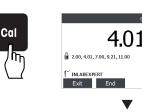

2013 14:34 /A 25.0°C MTC ©/Ā Value stable Cal 2

Press END to end calibration or repeat the last steps for a multi-point calibration.




Press SAVE to save calibration.



# SevenCompact<sup>™</sup> S210

### Switch display view

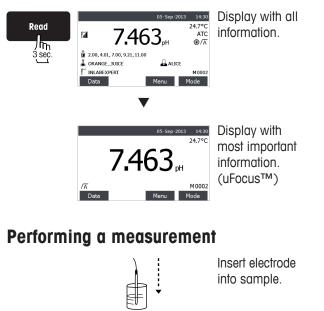



#### Mettler-Toledo AG, Analytical

CH-8603 Schwerzenbach, Switzerland Tel. +41 (0)44 806 77 11 Fax +41 (0)44 806 73 50 www.mt.com/pH

Subject to technical changes. © Mettler-Toledo AG 02/2014 30130406A



∕AValue stable

Endpoint type can be set to automatic, manual or timed in the menu settings.

#### Storing measurement data



Storage number is incremented.

Note: automatic storage to memory can be switched on in the setup menu.

#### Accessing menu settings



Press to enter the menu settings, e.g. to enter user ID, sample ID and Sensor ID.

