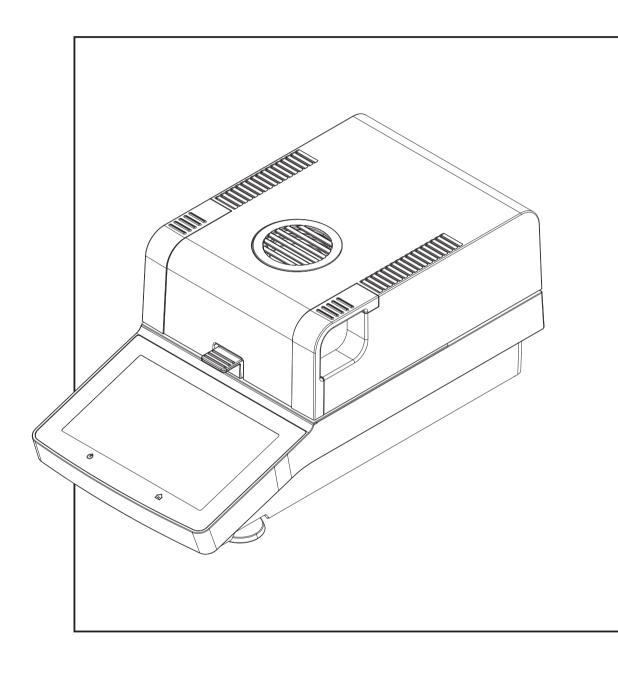
Moisture Analyzer

HC103



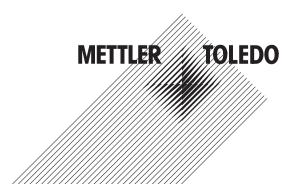


Table of Contents

1	Introd	uction	į
-	1.1	Further documents and information	Ę
	1.2	Explanation of conventions and symbols used	į
	1.3	Compliance information	(
2	Safety	Information	_
-	2.1	Definition of signal words and warnings symbols	-
	2.2	Product-specific safety notes	-
3	_	n and Function	10
	3.1	Overview	10
		3.1.1 Overview	10
		3.1.1.1 Drying unit	10
	2.0	3.1.2 Overview type plate	1
	3.2	Function description	1.
4	Install	lation and Putting into Operation	12
	4.1	Selecting the location	12
	4.2	Unpacking	12
	4.3	Scope of delivery	12
	4.4	Connecting the instrument	13
	4.5	Setup wizard	13
	4.6	Setting up the instrument	14
	4.7	Leveling the instrument	14
	4.8	Anti-theft device	15
	4.9	Setting date and time	16
	4.10	Adjusting after setup	16
	4.11	Transporting, packing and storing	16
		4.11.1 Transporting the Moisture Analyzer	16
			17
	110	4.11.3 Packing and storing	17
	4.12	Installing OSD device driver	
5		ting Elements and Interfaces	18
	5.1	Overview operation keys	18
	5.2	Electrical interfaces	18
6	Opera	tion	20
	6.1	Operation orientation	20
	6.2	User interface	20
		6.2.1 Operation keys	20
		6.2.2 Screen navigation	20
		6.2.3 Basic elements on the touch screen	2
		6.2.4 Icons and buttons	2
		6.2.4.1 Title bar icons	2
		6.2.4.2 Action bar buttons	2
		6.2.4.3 Status message icons	2
		6.2.5 Home screen	22
		6.2.6 Work screen	23
		6.2.7 Input dialogs	23
		6.2.7.1 Entering characters and numbers	23
		6.2.7.2 Entering numerical values	24
		6.2.7.3 Changing date and time	25
	6.3	6.2.8 Lists and tables	25
	C C	r coontinuo o altitule tricoatrement	_ / r

7	Softw	are Descri	ption	30
	7.1	Method de	efinition	30
		7.1.1	Method parameters	30
		7.1.1.1	Main measurement parameters	30
		7.1.1.2	Result and value handling	36
		7.1.1.3	Workflow handling	37
		7.1.1.4	General method properties	37
		7.1.2	Creating a method	38
		7.1.2.1	Creating a method with the Method Wizard	38
		7.1.2.2	Creating a method manually	40
		7.1.2.3	Creating a method by consulting the method library	40
		7.1.3	Editing a method	40
	7.2	Measuren	nent	41
		7.2.1	Performing a measurement	41
		7.2.2	Optimum sample preparation	44
		7.2.3	Work with weighing-in aid	44
		7.2.4	Working with shortcuts	44
	7.3	Results		45
		7.3.1	Graphical evaluations of measurement results	45
		7.3.2	Method versioning	48
		7.3.3	Exporting results	48
		7.3.3.1	Exporting results to a storage device	48
		7.3.3.2	Exporting results with EasyDirect Moisture	49
		7.3.3.3	Export settings	49
	7.4		st	50
		7.4.1	Tests	51
		7.4.1.1	Temperature test	51
		7.4.1.2	Weight test - external	53
		7.4.1.3	SmartCal test	53
		7.4.1.4	Measures after failed test	55
		7.4.2	Adjustments	55
		7.4.2.1	Weight adjustment - external	55
		7.4.2.2	Temperature adjustment	56
	7.5	7.4.3	History	58
	7.5	-	Hear profession	59
		7.5.1	User preferences	59
		7.5.1.1 7.5.1.2	Languages	59 60
		7.5.1.2	Screen	
		7.5.1.3	Sound	60 60
		7.5.1.4	Protection	61
		7.5.2.1	Instrument settings	61
		7.5.2.1	Peripherals	61
		7.5.2.2	Touch screen adjustment	63
		7.5.2.3	Applications settings	63
		7.5.3.1	Identifications management	63
		7.5.3.2	Print and export management	64
		7.5.4	Quality management	66
		7.5.4.1	Test / adjustment settings	66
		7.5.5	System and data management	70
		7.5.5.1	Export / Import	70
		7.5.5.2	Back up / Restore	71
		7.5.5.3	Export event logs	72
		7.5.5.4	Reset	72
		7.5.5.5	Update	72
		7.5.6	Help and tutorials.	73

8	Maint	enance	74
	8.1	Maintenance tasks	74
	8.2	Cleaning	74
		8.2.1 Sample chamber	75
		8.2.2 Heating module	
		8.2.3 Fan grill	
		8.2.4 Putting into operation after cleaning	
	8.3	Replacing the dust filter	77
	8.4	Replacing the power line fuse	77
	8.5	Updating the software	77
9	Troub	leshooting	79
	9.1	Error messages	79
	9.2	Error symptoms	79
10	Techn	ical Data	82
	10.1	General data	82
	10.2	Model-specific data	83
	10.3	Dimensions	84
	10.4	Interfaces specification	85
		10.4.1 RS232C	85
		10.4.2 USB host	85
		10.4.3 USB device	86
11	Acces	sories and Spare Parts	87
	11.1	Accessories	87
	11.2	Spare parts	90
12	Dispo	sal	92
	Index		93

3

Table of Contents

Moisture Analyzer

1 Introduction

Thank you for purchasing a Halogen Moisture Analyzer from METTLER TOLEDO.

Your Moisture Analyzer is fast, reliable and offers a high level of operating convenience and useful functions to facilitate determination of the moisture content of your samples.

The Moisture Analyzer conforms with all common standards and directives. It supports stipulations, work techniques and result records as demanded by all international quality assurance systems, e.g. GLP (Good Laboratory Practice), GMP (Good Manufacturing Practice). The instrument has a CE declaration of conformity and METTLER TOLEDO as the manufacturer has been awarded ISO 9001 and ISO 14001 certification. This provides you with the assurance that your capital investment is protected in the long term by a high product quality and a comprehensive service package (repairs, maintenance, servicing, adjustment service).

This document is based on the software version V 2.10.

EULA

The software in this product is licensed under the METTLER TOLEDO End User License Agreement (EULA) for Software.

www.mt.com/EULA

When using this product you agree to the terms of the $\ensuremath{\mathsf{EULA}}.$

1.1 Further documents and information

www.mt.com/moisture

This document is available in other languages online.

www.mt.com/HC103-RM

Instructions for cleaning a balance: "8 Steps to a Clean Balance"

www.mt.com/lab-cleaning-guide

Search for moisture software

www.mt.com/moisture-software

Search for documents

www.mt.com/library

For further questions, please contact your authorized METTLER TOLEDO dealer or service representative.

www.mt.com/contact

1.2 Explanation of conventions and symbols used

Conventions and symbols

Key and button designations are indicated by a picture or bold text (e.g. **Save**).

■ Note

For useful information about the product.



Refers to an external document.

Elements of instructions

In this manual, step-by-step instructions are presented as follows. The action steps are numbered and can contain prerequisites, intermediate results and results, as shown in the example. Sequences with less than two steps are not numbered.

- Prerequisites that must be fulfilled before the individual steps can be executed.
- 1 Step 1
 - Intermediate result
- 2 Step 2
- → Result

Moisture Analyzer Introduction

1.3 Compliance information

European Union

The instrument complies with the directives and standards listed on the EU Declaration of Conformity.

United States of America

This equipment has been tested and found to comply with the limits for a **Class A** digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

The FCC Supplier Declaration of Conformity is available online.

► http://www.mt.com/ComplianceSearch

Introduction Moisture Analyzer

2 Safety Information

Two documents named "User Manual" and "Reference Manual" are available for this instrument.

- The User Manual is printed and delivered with the instrument.
- The electronic Reference Manual contains a full description of the instrument and its use.
- Keep both documents for future reference.
- Include both documents if you transfer the instrument to other parties.

Only use the instrument according to the User Manual and the Reference Manual. If you do not use the instrument according to these documents or if the instrument is modified, the safety of the instrument may be impaired and Mettler-Toledo GmbH assumes no liability.

2.1 Definition of signal words and warnings symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

Signal words

DANGER A hazardous situation with high risk, resulting in death or severe injury if not avoided.

WARNING A hazardous situation with medium risk, possibly resulting in death or severe injury if

not avoided.

CAUTION A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.

NOTICE A hazardous situation with low risk, resulting in damage to the instrument, other

material damage, malfunctions and erroneous results, or loss of data.

Warning symbols



General hazard: read the User Manual or the Reference Manual for information about the hazards and the resulting measures.



Hot surface



Notice

7

2.2 Product-specific safety notes

Intended use

This instrument is designed to be used by trained staff. The instrument is intended for determining the weight loss during drying of samples.

Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

Moisture determination applications must be optimized and validated by the user according to local regulations. Application-specific data provided by METTLER TOLEDO is intended for guidance only.

Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties.

Mettler-Toledo GmbH assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. Mettler-Toledo GmbH assumes that the instrument owner provides the necessary protective gear.

Moisture Analyzer Safety Information

Personal protective equipment



Insulated gloves



Lab coat



Goggles

Safety notes



↑ WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to injury and death. If the instrument cannot be shut down in an emergency situation, people can be injured and the instrument can be damaged.

- 1 Check that the voltage printed on the instrument is the same as your local power supply voltage. If this is not the case, under no circumstances connect the instrument to the power supply, but contact a METTLER TOLEDO representative.
- 2 Only use the three-core power cable with equipment grounding conductor supplied by METTLER TOLEDO to connect your instrument.
- 3 Only connect it to a three-pin power socket with earthing contact.
- 4 Only use standardized extension cables with equipment grounding conductor for operating the instrument.
- 5 Make sure that the power plug is accessible at all times.
- 6 Arrange the cables so that they cannot be damaged or interfere with the operation.
- 7 Keep all electrical cables and connections away from liquids.



⚠ WARNING

Injury or death due to toxic or corrosive substances

Heating up toxic or corrosive substances, e.g., acids, can result in toxic or corrosive vapors that can cause injuries if they come in contact with the skin or the eyes or if they are inhaled.

- 1 When using chemicals and solvents, comply with the instructions of the manufacturer and the general laboratory safety rules.
- 2 Set up the instrument in a well-ventilated location.
- 3 When using dry substances that form toxic gases, place the instrument in a fume hood.



WARNING

Death or serious injuries due to flammable solvents

Flammable solvents in the vicinity of the instrument can ignite and lead to fire and explosions.

- 1 Keep flammable solvents away from the instrument.
- When using chemicals and solvents, comply with the instructions of the manufacturer and the general laboratory safety rules.

Safety Information Moisture Analyzer

555

CAUTION

Burns due to hot surfaces

During operation, parts of the instrument can reach temperatures that can cause burns if touched.

- 1 Do not touch the area marked with the warning symbol.
- 2 Ensure sufficient free space around the instrument to avoid heat accumulation and overheating (approx. 1 m free space above the heating module).
- 3 Never cover, tape or clog the vent over the sample chamber. Do not tamper with the vent in any other way.
- 4 Exercise caution when removing a sample. The sample itself, the sample chamber, the draft shield and the sample pan may be very hot.
- 5 Do not open the heating module during operation. Always let it cool down completely before opening.
- 6 Do not modify the heating module in any way.



NOTICE

Damage to the instrument due to corrosive substances and vapors

Corrosive substances and corrosive vapors can damage the instrument.

- 1 When using chemicals and solvents, comply with the instructions of the manufacturer and the general laboratory safety rules.
- 2 Ensure that the instrument parts touching your sample substance can not get altered by it.
- 3 Wipe off any condensation of corrosive vapours after an operation.
- 4 Work with small samples.



NOTICE

Damage to the instrument or malfunction due to the use of unsuitable parts

 Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

Moisture Analyzer Safety Information


9

3 Design and Function

3.1 Overview

3.1.1 Overview

3.1.1.1 Drying unit

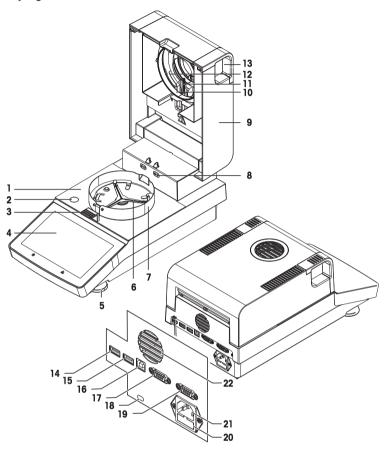

Legend

- 1 Sample chamber
- 3 Sample pan handler
- 5 Leveling screw
- 7 Draft shield element
- **9** Heating module (with thermal overload protection)
- **11** Temperature sensor
- 13 Handles for opening the sample chamber
- **15** USB Host 2
- 17 RS232C

10

- **19** RS232C (for production only)
- 21 Power supply socket


- 2 Level indicator
- 4 Touch screen
- 6 Sample pan holder
- 8 Contacts for temperature adjustment kit (optional)
- 10 Reflector ring
- 12 Protective glass
- **14** USB Host 1
- 16 USB Device
- **18** Kensington slot for anti-theft purposes
- 20 Power line fuse
- **22** Fan

Design and Function Moisture Analyzer

3.1.2 Overview type plate

Drying unit type plate

The type plate of the drying unit is located on the side of the drying unit and contains the following information:

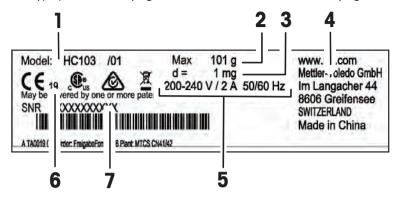


- 1 Model designation
- 3 Maximum capacity
- **5** Power supply
- **7** Serial number (SNR)

- 2 Readability
- 4 Manufacturer
- 6 Year of manufacture

3.2 Function description

METTLER TOLEDO Moisture Analyzers work on the thermogravimetric principle. At the start of the measurement the Moisture Analyzer determines the weight of the sample, the sample is then quickly heated by the absorption of the emitted infrared radiation. During the drying process the instrument continually measures the weight of the sample and displays the weight reduction, up to the final result.

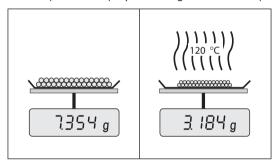

With halogen heating technology, maximum heating power is quickly reached and allows use of high temperatures. Uniform heating of the sample material ensures repeatability of the drying results and makes it possible to use small amounts of sample.

The METTLER TOLEDO Moisture portfolio comprises a range of moisture analyzers that differ from each other in both hardware and software.

The following features are available for the HC103 model:

- Robust and compact design, ideal for limited laboratory space.
- Intuitive operation thanks to a large touchscreen and easy to navigate user interface. See [User interface ▶ Page 20]
- Quick and easy cleaning. See [Cleaning ▶ Page 74]
- Quick access to methods and results with shortcuts on the home screen. See [Working with shortcuts > Page 44]
- Easy reporting. See [Exporting results ▶ Page 48]
- ID management. See [Identifications management ▶ Page 63]
- Setup wizard and instrument tutorial. See [Setup wizard ▶ Page 13] and [Help and tutorials ▶ Page 73]

11

Moisture Analyzer Design and Function

4 Installation and Putting into Operation

4.1 Selecting the location

Requirements of the location

Place indoors on stable table

Ensure sufficient spacing Level the instrument provide adequate lighting table

Avoid direct sunlight

Avoid vibrations

Avoid strong drafts

Avoid temperature fluctuations

Sufficient distance for Moisture Analyzers: > 15 cm next to the instrument, > 1 m above the heating module.

4.2 Unpacking



Note

Store all parts of the packaging. This packaging guarantees the best possible protection for the transport of your instrument.

After receiving the instrument, proceed as follows:

- 1 Open the package and remove drying unit and the accessories.
- 2 Remove the packaging from the instrument.
- 3 Check the Moisture Analyzer for transport damage and complete scope of delivery.
- 4 In the event of damage or missing accessories, immediately inform a METTLER TOLEDO representative.

4.3 Scope of delivery

Moisture Analyzer	Documentation	Accessories
1 Drying unit with terminal	1 User Manual	80 aluminium sample pans
• 1 Draft shield	1 Application brochure «Guide	3 Specimen samples (glass)
• 1 Power cable	to Moisture Analysis»	fibre filter)
1 Sample handler	EU-countries: 1 CE declaration	1 SmartCal sampler
1 Sample pan holder	of conformity	

4.4 Connecting the instrument

Connecting the instrument



⚠ WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to injury and death. If the instrument cannot be shut down in an emergency situation, people can be injured and the instrument can be damaged.

- 1 Check that the voltage printed on the instrument is the same as your local power supply voltage. If this is not the case, under no circumstances connect the instrument to the power supply, but contact a METTLER TOLEDO representative.
- 2 Only use the three-core power cable with equipment grounding conductor supplied by METTLER TOLEDO to connect your instrument.
- 3 Only connect it to a three-pin power socket with earthing contact.
- 4 Only use standardized extension cables with equipment grounding conductor for operating the instrument.
- 5 Make sure that the power plug is accessible at all times.
- 6 Arrange the cables so that they cannot be damaged or interfere with the operation.
- 7 Keep all electrical cables and connections away from liquids.

Two different versions of drying units with country-specific power cable are available:

110 V AC or 230 V AC

- Instrument is at the final location.
- 1 Connect the power cable to the power supply socket on the instrument.
- 2 Connect the power cable to the power supply.

See also

Drying unit ▶ Page 10

4.5 Setup wizard

When you switch on your Moisture Analyzer for the first time or when a factory reset was performed, the **Setup Wizard** appears. The **Setup Wizard** is a step-by-step instruction, that helps to define the basic functions (**Language** and **Date format**) and to install the sample chamber of your device. The buttons in the action bar can be used for navigation.

It must be successfully finished. If it is not finished, the device is not operational and the function will reappear when the device is switched on the next time.

After the application **Setup Wizard** is successfully finished, the settings can be changed via **Settings** of the user home screen.






Note


The application **Instrument tutorial** starts automatically after the application **Setup Wizard** has successfully finished. This option can be skipped by inactivating the checkbox **Proceed to the Instrument Tutorial**. If you want to redo the application **Instrument tutorial**, it can be found in the section **Settings**.

4.6 Setting up the instrument

- Instrument is connected to the power supply.
- 1 Open the sample chamber.
- 2 Place the draft shield element. Place the notches onto the screw heads on the bottom of the sample chamber.

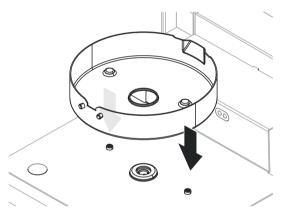

- 3 Carefully insert the sample pan holder. Make sure that the sample pan holder is correctly positioned and falls into position (see picture).
- 4 Insert the sample pan handler.
- 5 Press [6] to start the instrument.

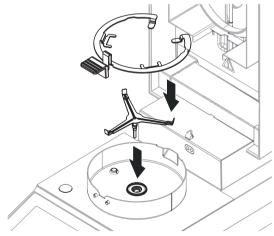

4.7 Leveling the instrument

Exact horizontal positioning and stable installation are prerequisites for repeatable and accurate results. To compensate for small irregularities or inclinations $(\pm 2\%)$ at the location, the instrument must be leveled.



Note

For leveling the drying unit use the function **Leveling guide** that can be found on **Settings** > **Help and tutorials** > **Instrument tutorial** > **1. Leveling the instrument**.

For exact horizontal positioning, the device has a level indicator and two leveling feet. When the air bubble in the level indicator is exactly in the center, the instrument is standing perfectly horizontally.



Note


The drying unit should be re-leveled each time its location is changed.

To level it, proceed as follows:





- 1 Position your Moisture Analyzer at the selected location.
- 2 Turn the two leveling feet until the air bubble is in the center of the level indicator.

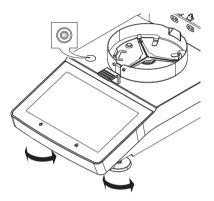

Air bubble at 12 o'clock:



turn both feet clockwise.



Air bubble at 3 o'clock:



turn left foot clockwise, right foot counterclockwise.



Air bubble at 6 o'clock:



turn both feet counterclockwise.




Air bubble at 9 o'clock:



turn left foot counterclockwise, right foot clockwise.

4.8 Anti-theft device

For protection against theft, the Moisture Analyzer is fitted with a Kensington security slot for an anti-theft device. To find a suitable anti-theft cable **see** [Accessories Page 87].

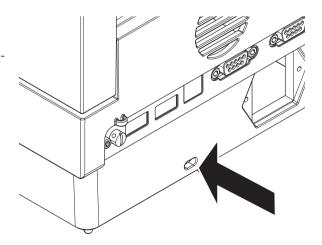

4.9 Setting date and time

Navigation: Home > Settings > Instrument settings > Regional Settings

When the instrument was switched into operation for the first time, date, time and language have been defined with the function **Setup Wizard**. These settings are retained even if the instrument is being disconnected from the power supply. The settings can also be changed manually as follows:

Set the current date

- Regional Settings is selected.
- 1 Tap Date.
- 2 Set day, month and year.
- 3 Confirm with Set date.

Set the current time

- Regional Settings is selected.
- 1 Tap **Time**.
- 2 Set hours and minutes.
- 3 Confirm with Set time.

4.10 Adjusting after setup

To obtain accurate measuring results, it is necessary to adjust the integrated balance as well as the heating module under working condition.

Adjusting is necessary before the instrument is used for the first time or after a change of location.

The following adjustment options are possible:

- Balance adjustment with external weight
- Temperature adjustment with temperature kit
- To verify the overall performance of the Moisture Analyzer after the aforementioned adjustments, perform a SmartCal test.


For more information to perform these operations **see** [Test/Adjust ▶ Page 50].

4.11 Transporting, packing and storing

4.11.1 Transporting the Moisture Analyzer

Transporting the Moisture Analyzer over short distances

- 1 Switch off the instrument and unplug all interface cables.
- 2 Hold the instrument with both hands as shown.
- 3 Carefully lift the instrument and carry it in horizontal position to the new location.

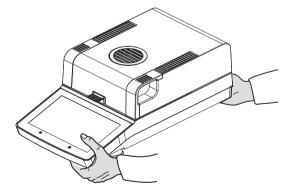

Transporting the Moisture Analyzer over long distances

For transporting the Moisture Analyzer over long distances, always use the original packaging or the designated transport case (**see** "Accessories").

4.11.2 Putting into operation after transport

Putting into operation after transport:

- 1 Connect the instrument to the power supply.
- 2 Check the level status. Level the Moisture Analyzer if necessary.
- 3 METTLER TOLEDO recommends to perform weight and temperature tests and, if necessary, adjustments after transporting the Moisture Analyzer.
- The Moisture Analyzer has been put into operation and is ready to use.

See also

- Connecting the instrument ▶ Page 13
- Test/Adjust ▶ Page 50

4.11.3 Packing and storing

Packaging

Store all parts of packaging in a safe place. The elements of the original packaging are developed specifically for the Moisture Analyzer and its components to ensure optimal protection during transportation or storing.

Storing

Store the Moisture Analyzer under following conditions:

- Indoor and in the original packaging.
- According to the environmental condition, see "Technical data".

Note

When storing for longer than six months, the rechargeable battery may be down (date and time get lost).

4.12 Installing USB device driver

To transmit commands via USB device interface, a USB driver program has to be installed on the instrument and PC. The driver program is automatically installed during setup of the instrument. However, in case of an installation error, the software can be downloaded online and installed manually.

Find the download link on:

www.mt.com/labweighing-software-download

Requirements

- PC with one of the following Microsoft Windows® 32-bit/64-bit operating systems: Win 7 (SP1), Win 8 or Win 10.
- USB cable to connect the instrument to the PC.

Downloading and installing the USB driver


- 1 Open the download portal (see above) on your PC.
- 2 Download the driver software and extract the .zip file.
- 3 Right-click on the extracted .exe installation program and select **Run as Administrator**.
- 4 If a safety warning appears, allow windows to perform the installation.
- 5 Click **Next** and follow the installer's instructions.

Connecting the instrument to the PC

- 1 Connect the moisture analyzer to the preferred USB port on the PC.
- 2 Follow the instructions on the PC and install the software automatically (recommended).
- 3 Check that the instrument is listed as a device (COM Port) on your PC.
- 4 Set USB interface in the peripheral settings of the instrument.


5 Operating Elements and Interfaces

5.1 Overview operation keys

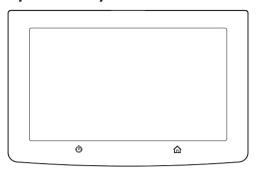

Ф	û
ON/OFF	Home

5.2 Electrical interfaces

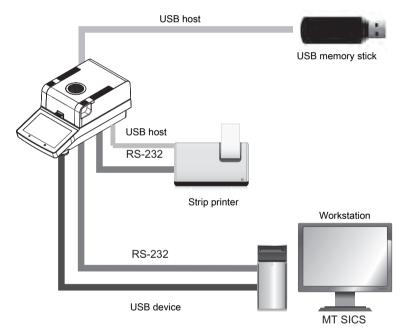

Note

The USB host connection can only be used with specific devices. The following devices are officially supported:

- USB memory stick (for specific characteristics see list below)
- A bar code reader can be connected to the USB port. With a connected bar code reader, bar codes can be added in a keyboard dialog.

Note for using USB memory stick

- USB memory sticks up to 32 GB are supported.
- Recommended formatting: FAT32 (NTFS not supported).
- Max. cluster size: 32 KB.
- Max. file size: 32 MB.

- Make sure that the stick is fully inserted.
- Make sure that the write protection is deactivated if present.

6 Operation

6.1 Operation orientation

The Moisture Analyzer is mainly operated via the touchscreen and keys of the terminal.

Before a measurement can be made, a method must first be defined. Methods define the parameters of a measurement, such as the temperature to which the sample is heated or the criterion after which the measurement is finalized. The optimum parameters differ from sample to sample. The instrument and METTLER TOLEDO support you in finding a suitable method for your sample.

Once a method has been defined, the actual measurement can be started: The sample is placed in the sample chamber and prepared. Then the heating module is closed and the sample is heated. Depending on the sample and the method, the measurement can take from a few minutes to a few hours.

When the measurement is finished, the end result is shown on the terminal screen. The result of the measurement can also be exported or printed.

The Moisture Analyzer can be used immediately for another measurement.

6.2 User interface

6.2.1 Operation keys

K	Yey	Designation	Description
	ψ		Switches the Moisture Analyzer on / into standby mode (to switch it off completely, it has to be unplugged from the power supply).
		Home	Directly returns to the home screen from any other menu level.

6.2.2 Screen navigation

The colored touch screen is a touch-sensitive WVGA monitor. It displays information and allows you to enter commands by touching certain areas on its surface: You can choose the information displayed on the screen, change terminal settings or perform certain operations on the instrument.

Only those elements which are available for the current dialog appear on the display.



NOTICE

Damage to the touch screen due to pointed or sharp objects

Operate the touch screen with your fingers.

Note

The sensitivity of the touch screen is set to a default value. The sensitivity of the touch screen can be defined with the function **Touch screen adjustment** in the section **Settings**.

The surface navigation of the touch screen works in the same way as on most common touch screens:

Selecting a button or an icon

- Tap on it.

Scrolling

- Drag the button in the scroll bar up or down.
- Or tap ▲ or ▼

Operation Moisture Analyzer

6.2.3 Basic elements on the touch screen

The graphical user interface consists of the following basic elements.



	Name	Explanation
1	Status message icons	Show information about the instrument status. The quick access icon (for settings of brightness and sound) is displayed continuously. See Status Message Icons
2	Status bar	Shows the user name, status icons as well as date and time.
3	Title bar	Shows title of the current screen, contains special functions. See [Title bar icons ▶ Page 21].
4	Content area	Main working area of the menus and the applications. Shows content depending on application or performed action. May also show graphic illustrations such as the curve diagram of a data set (e.g. drying curve).
5	Action bar	Contains action buttons which are required and available in the current dialog. See [Action bar buttons ▶ Page 21].

6.2.4 Icons and buttons

6.2.4.1 Title bar icons

The title bar icons only appear when they are active. The icons which appear in the title bar, are listed with their function below:

Parameter	Description
	Create a shortcut and add it to your user home screen. Edit or remove an existing shortcut see Using Shortcuts.

6.2.4.2 Action bar buttons

The action bar contains action buttons for triggering any of the actions which are required and available in the current dialog (e.g. **Back**, **->0/T<-**, **Print**, **Save**, **Delete**, **OK**). There are up to six action buttons available for a work step.

6.2.4.3 Status message icons

Status messages are displayed by means of small icons in the status bar of the display. The status icons indicate the following:

Icon	Status description	Diagnostic	Remedy
Se Se	Service is due.	_	Contact your METTLER TOLEDO support representative.

21

Moisture Analyzer Operation

Icon	1	Status description	Diagnostic	Remedy
)	ď	Quick Access icon (displayed continuously)	Quick setting of brightness and sound.	-
		EasyDirect Moisture connection icon	Indicates the connection of the device to EasyDirect Moisture .	-

6.2.5 Home screen

22


The user **Home** screen is the main screen and appears after startup or login of the instrument. All screens of the user interface can be accessed from the home screen. Returning to the **Home** screen is possible from all other screens of the user interface by pressing the [$\upphantom{\up$



	Name	Explanation
1 Main Menu Measurement		Measurement
Start measurement.		Start measurement.
		Prerequisite: Method is defined and all necessary settings are made.
		Results
		Display, print and export results
		Method Definition
Define, edit, test or delete a method		Define, edit, test or delete a method
		Test/Adjust
		Adjust or test the integrated balance and the heating module and perform SmartCal tests.
		Settings
Define instrument-, user-, and data-management settings. Find help under this menu item.		Define instrument-, user-, and data-management settings. Find help and tutorials under this menu item.
2	User shortcuts	Shows user-specific shortcuts for frequently used methods. Shortcuts are saved in the user profile.
3	Instrument info	Shows general info about the instrument and the software (e.g. serial number, software version).

Operation Moisture Analyzer

6.2.6 Work screen

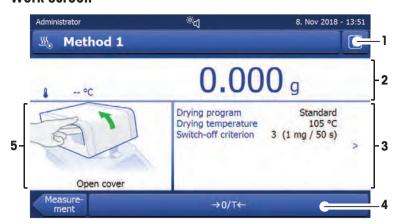

	Name	Explanation	
1	Shortcut button	Adding/editing a shortcut to the current method (to the home screen).	
2	Value panel	Current measured (or predicted) values of the work process.	
3	Parameter panel	Parameters of the current work process. A detailed overview of the method parameters appears after tapping the parameter panel.	
		Identifications (IDs) appear after tapping the ID panel for entering or editing values (comments). ID panel appears only if identification input is activated in the menu.	
4	Action buttons	Action buttons which are required and available for the current dialog (e.g. Back , -> 0/T<- , Print , Save , Delete , OK).	
5	Graphics panel	Graphical illustrations, e.g. of the drying curves, instructions for the user for performing tasks, and weighing-in aid.	

6.2.7 Input dialogs

6.2.7.1 Entering characters and numbers

The keyboard dialog allows you to enter characters including letters, numbers and a variety of special characters. The country-specific layout of the keyboard is in accordance with the selected language, **see** Regional Settings.



	Name	Explanation
1	Input field	Shows the data that has been entered.
2	Explanation field	Indicates the maximum number of characters allowed.
3	Shift	Switches the keyboard between standard and special characters.

23

Moisture Analyzer Operation

- Confirm with **[OK]** (To exit keyboard tap **[Cancel]**).

	Function
슢	Shifts between lower and upper case letters
\boxtimes	Deletes last character
←	Cursor left
\rightarrow	Cursor right

Note

It is also possible to place cursor directly in input field by tapping the respective position.

6.2.7.2 Entering numerical values

The numerical keypad dialog allows you to enter numerical values.



	Name	Explanation
1	Switch	If a switch appears, located on the left side in the title bar; it must be switched on for the entry of a value: $O = \mathbf{OFF}$, $I = \mathbf{ON}$
2	Input field	Shows the data that has been entered.
3	Explanation field	The valid input range is indicated.
\boxtimes	Delete button	Deletes last character.
(Cursor leff	Jumps one character to the left.
\rightarrow	Cursor right	Jumps one character to the right.


- Confirm with [**OK**].

To exit keypad, tap [Cancel].

Operation Moisture Analyzer

6.2.7.3 Changing date and time

This dialog (Picker view) allows you to set the date and time by picking from the built-in calendar/clock.

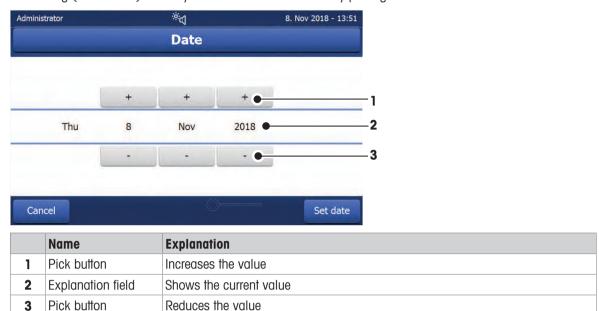

- Confirm with [Set date]. (To exit without changing, tap [Cancel].)

6.2.8 Lists and tables

Simple list view

Basic elements of a simple list including a content title. If necessary, the content area can be scrolled vertically.



	Name	Explanation	
1	Content title	Shows the subject of the current content	
2	Content area	Shows the relevant content	
3	Arrow icon	[♣] Shows the list in ascending order.	
		[1] Shows the list in descending order.	
4	Scrollbar	oar [🛕] Scroll up	
5		Scroller	
6		[▼] Scroll down	

Moisture Analyzer Operation

25

Further types of lists

Accordion panels are opened or closed by tapping [\alpha].



Closable panels are opened or closed by tapping [▶].



6.3 Performing a simple measurement

To check the function of the instrument and to become familiar with the process, METTLER TOLEDO recommends carrying out a simple example measurement, e.g. using a glass fibre filter to measure the moisture content of water. During this first measurement, the instrument works with factory default settings.

- The instrument is connected to the power supply.
- To switch on, press 心.
- 1 Tap **Method Definition**.
 - → Method Definition menu appears.
- 2 To define a new method, tap **New...**.
- 3 To define a new method manually, tap **Define method manually**.
 - → The keypad appears.
- 4 Enter a name for your first method e.g. **My first method**.
- 5 Confirm with **OK**.
- 6 To store your new method with factory settings, tap **Save**.
- 7 Tap Home.
- User home screen appears.



26

Note

The number of methods that can be created is limited to 20.

Operation Moisture Analyzer


Select the method for measuring

- 1 Tap **Measurement**.
 - → The methods list is displayed.
- 2 Tap My first method.
 - → The work screen of the method My first method appears.
- 3 Open the sample chamber.

Placing the accessories

- The display now prompts you to load the predefined accessories and tare the balance.
- 1 Place the empty accessories in the sample pan handler.
- 2 Place the sample pan handler in the sample chamber. Ensure that the tongue of the pan handler fits exactly in the slot of the draft shield element. The sample pan must lie flat in the pan holder.

We advise you to work with the sample pan handler at all times. The pan handler is ergonomic, automatic positioning, safe and provides protection against possible burns due to the hot accessories.

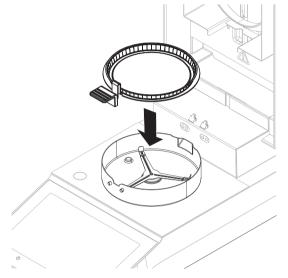

27

Taring the balance



Note

Before the sample pan is placed into the sample chamber, the left corner of the display must show **Open cover**, place sample pan and tare.

If the display shows **Add sample**, tap **Cancel** before placing the empty sample pan in the sample chamber.


- 1 Close the sample chamber.
 - → The instrument tares the balance (Starting mode: Automatic).
- 2 After tarring, open the sample chamber.

Note

Moisture Analyzer Operation


In this example, the glass fiber filter is part of the sample for the sake of simplicity. In routine liquid measurements, the glass fiber filter is not part of the sample and must be tared together with the sample pan.

1 Place the specimen sample in the sample pan.

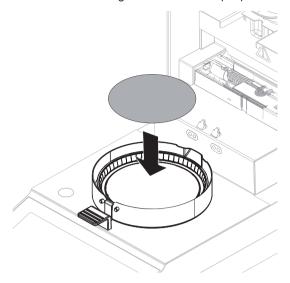

- 2 Wet the specimen sample with a few drops of water.
- 3 Close the sample chamber.
 - → The drying process starts automatically.

Drying process

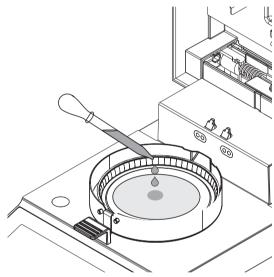

You can follow the measurement process on the display.

- The drying process is continuously displayed graphically.
- The current temperature in the heating module is displayed as well as the elapsed drying time and the current drying value.
- The display shows the selected settings.
- Tap on Stop drying. The options Abort measurement and save data or Cancel are available.
- Tap on Abort measurement and save data to abort the process.
- Tap on Cancel to continue the process.

At the end of the drying process you can read the moisture content of your sample on the display.



CAUTION

Burns due to hot surfaces!

Parts of the instrument can reach temperatures that cause injuries if touched. Sample, sample pan and other parts inside the sample chamber may still be hot.

Do not touch the place marked with the warning symbol.


Operation Moisture Analyzer

- The drying process is completed.
- 1 Open the sample chamber.
- 2 Carefully remove the sample pan handler from the sample chamber.

Note

To remove the sample pan from the handler, lift the pan slightly from below and pull it out of the handler.

- 3 To perform a further measurement with the current method, tap **Next Sample**.
 - To perform a measurement with a new method, tap **Methods**.
 - To return to the home screen, press 🔔.

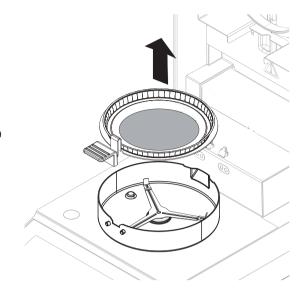

29

Moisture Analyzer Operation

7 Software Description

7.1 Method definition

Navigation: Home > Method definition

This function offers you to define new methods and to change, delete, export or import existing methods depending on the user rights. It allows you to store up to 20 individual methods.

All parameters used in a drying method can be defined under this menu item.

Resources on methods and method definition:

www.mt.com/moisture-guide

www.mt.com/moisture-methods

7.1.1 Method parameters

7.1.1.1 Main measurement parameters

Navigation: Home > Method Definition > Method name > Main measurement parameters

These parameters are relevant for measuring and must be determined for all samples.

The following parameters can be defined:

Menu item	Explanation	Further Information
Defines the drying program which is best suite the specific sample. (including temperature ar switch-off criteria)		see [Drying program ▶ Page 30]
Display mode	Defines the type of value for displaying and printing.	see [Display mode ▶ Page 34]
Start weight	Defines a start weight so that the samples always have approximately the same weight.	see [Start weight ▶ Page 35]

7.1.1.1.1 Drying program

Navigation: Home > Method Definition > Method name > Main measurement parameters > Drying program

This function offers you different predefined drying programs for optimum matching of the drying characteristics to the sample used. Furthermore, the drying temperature and the switch-off criterion can be set.



Note

You will find more information about defining methods in the delivered application brochure "Guide to Moisture Analysis".

The following parameters can be defined:

Parameter	Description	Values
Drying program	Defines the drying program which is best suited for the specific sample.	Standard* I Rapid

^{*} Factory setting

Further information:

- Standard drying see [Standard drying ▶ Page 31]
- **Rapid** drying **see** [Rapid drying ▶ Page 31]

Software Description Moisture Analyzer

Types of drying programs

Standard drying



Standard drying

This drying program is set in the factory and is suitable for most samples. The sample is heated to the drying temperature.

Factory setting

Drying temperature = $105 \, ^{\circ}\text{C}$, Switch-off criterion $3 = 1 \, \text{mg} / 50 \, \text{s}$

The following parameters can be defined:

Parameter	Description	Values
Drying temperature	Defines the drying temperature.	40230 °C (105 °C *)
Switch-off criterion	Defines the criterion when the instrument should end the drying.	1 (1 mg / 10 s) 2 (1 mg / 20 s) 3 (1 mg / 50 s)* 4 (1 mg / 90 s) 5 (1 mg / 140 s) Timed Free (mg / s) Free (% / s)

^{*} Factory setting

Setting the drying temperature

Navigation: Method Definition > Name of method > Main measurement parameters > Drying temperature In this menu item you can set the final drying temperature.



Note

The admissible input range for the drying temperature is indicated.



- 1 Start of drying
- 2 Final temperature

Setting switch-off criterion

see [Switch-off criterion (SOC) ▶ Page 32]

Rapid drying



Rapid drying

This drying program is primarily suitable **for samples with a moisture content over 30%**. Following the start, the selected temperature is exceeded by 40% for 3 minutes (the maximum heating temperature is 230 °C) to compensate for the cooling due to vaporization and to accelerate the drying process. The drying temperature is then lowered to the set value and maintained.

31

Factory setting

Drying temperature = $105 \, ^{\circ}\text{C}$, Switch-off criterion $3 = 1 \, \text{mg} / 50 \, \text{s}$

Moisture Analyzer Software Description

The following parameters can be defined:

Parameter	Description	Values
Drying temperature	Defines the drying temperature.	40230 °C (105 °C *)
Switch-off criterion	Defines the criterion when the instrument should end the drying.	1 (1 mg / 10 s) 2 (1 mg / 20 s) 3 (1 mg / 50 s)* 4 (1 mg / 90 s) 5 (1 mg / 140 s) Timed Free (mg / s) Free (% / s)

^{*} Factory setting


Setting the drying temperature

Navigation: Method Definition > Name of method > Main measurement parameters > Drying temperature In this menu item you can set the final drying temperature.



Note

The admissible input range for the drying temperature is indicated.

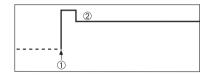

- 1 Start of drying
- 2 Final temperature

Setting switch-off criterion

see [Switch-off criterion (SOC) ▶ Page 32]

Switch-off criterion (SOC)

Navigation: Home > Method Definition > Method name > Main measurement parameters > Drying program

This function offers you different switch-off criteria. A switch-off criterion defines when the instrument should end the drying. Switch-off criteria ensure that measurements are finished in the same condition (weight loss per unit of time) every time, ensuring repeatable measurements.



Accurate results require same start weights, **see** [Start weight ▶ Page 35].

The following parameters can be defined:

Parameter	Description	Values
Switch-off criterion	Defines the criterion when the instrument should end the drying.	1 (1 mg / 10 s) 2 (1 mg / 20 s) 3 (1 mg / 50 s)* 4 (1 mg / 90 s) 5 (1 mg / 140 s) Timed Free (mg / s) Free (% / s)

^{*} Factory setting

- Weight loss per unit of time (5 preprogrammed settings)
- Free switch-off criterion (2 different settings)
- Timed switch-off

Software Description Moisture Analyzer

Weight loss per unit of time

The switch-off defines in which drying condition, the measurement is completed. This switch-off is based on a weight loss per unit of time. As soon as the mean weight loss is less that a preset value during a specified time, the instrument considers drying as complete and automatically discontinues the measurement process.




Note

The switch-off criterion is inactive during the first 30 sec.

The following 5 preprogrammed settings are available

Switch-off criterion	Δg	Δt	Description	
1	1 mg	10 s	This setting is suitable for fast measurements to determine a trend.	
2	1 mg	20 s	This setting is suitable for quick drying samples.	
3	1 mg	50 s	This is the factory setting . It is suitable for most kind of samples.	
4	1 mg	90 s	This setting is suitable for samples which dry moderately quickly or for samples with higher accuracy requirements.	
5	1 mg	140 s	This setting is suitable for samples which dry very slowly (trapped moisture, skin formation) or for samples with very low moisture content e.g. plastics. Not suitable for very temperature sensitive samples.	

The graph below exemplifies the mode of operation of the switch-off operation (not to scale).

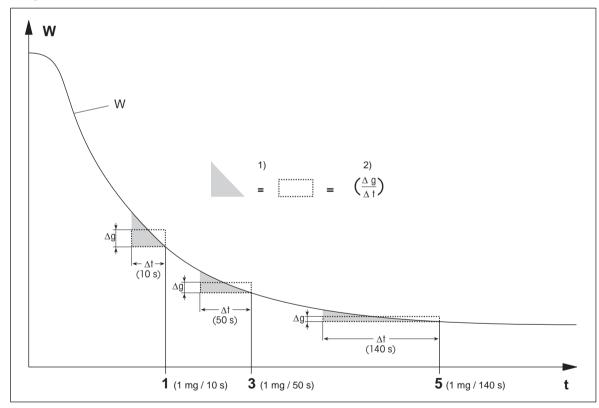

t = time

W

1, 3, 5

- = weight of sample
 - switch-off criteria shown as an example
- 1) = equal area
- 2) = mean weight loss per unit of time

33

Free switch-off criterion

The free switch-off criterion is based on a user-defined mean weight loss per unit of time. The following 2 settings are available:

- Switch-off criterion > Free (mg / s)... (weight loss per unit of time)
- Switch-off criterion > Free (% / s)... (weight loss in percent per unit of time)

Moisture Analyzer Software Description

Timed switch-off

With this switch-off criterion the measurement lasts until the pre-set drying time has elapsed. The display provides you with continuous information in the drying time.

7.1.1.1.2 Display mode

Navigation: Home > Method Definition > Method name > Main measurement parameters > Display mode

With this function you can select the desired type of result display. You also define what type of values are printed on the records.

The following parameters can be defined:

Parameter	Description	Values
Display mode	Defines the type of value for displaying and printing.	%MC* %DC %AM %AD g -%MC

^{*} Factory setting

Explanation of the values

- **%MC** Moisture Content (calculated value)
- %DC Dry Content (calculated value)
- **%AM** ATRO Moisture Content (calculated value)
- %AD ATRO Dry Content (Wet weight, calculated value)
- **g** Weight in grams
- -%MC Moisture Content (negative value, calculated value)



Note


Calculated values are indicated with an asterisk in the display.

Detailed information:

%MC - Moisture Content

The moisture content of the sample is displayed (and printed out) as a percentage of the wet weight ($WW = initial\ weight = 100\%$). This is the **factory setting**.

During the measurement the value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "%MC" (Moisture Content, e.g. 11.35 %MC) also for the printed results.

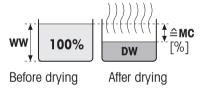

$$MC = \frac{WW - DW}{WW} \cdot 100 \%$$

MC = Moisture Content [0...100%]


WW = wet weight

DW = dry weight

%DC - Dry Content

The dry content of the sample is displayed (and printed out) as a percentage of the wet weight (WW = initial weight = 100%).

During the measurement the value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "%DC" (Dry Content, e.g. 88.65 %DC) also for the printed results.

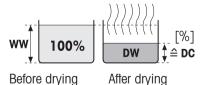

$$DC = \frac{DW}{WW} \cdot 100 \%$$

DC = dry content [100...0%]


WW = wet weight

DW = dry weight

%AM - ATRO Moisture Content 1)

The moisture content of the sample is displayed (and printed out) as a percentage of the dry weight (DW = final weight = 100%)

During the measurement the value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "%AM" (ATRO Moisture Content, e.g. 255.33 %AM) also for the printed results.

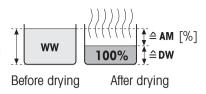

$$AM = \frac{WW - DW}{DW} \cdot 100 \%$$

AM = ATRO moisture content [0...1000%]

WW = wet weight

DW = dry weight

%AD - ATRO Dry Content (Wet weight)1)

The wet weight of the sample is displayed (and printed out) as a percentage of the dry weight (DW = final weight = 100%)

During the measurement the value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "%AD" (ATRO Dry Content, e.g. 312.56 %AD) also for the printed results..

$$AD = \frac{WW}{DW} \cdot 100 \%$$

AD = ATRO dry content [100...1000%]

WW = wet weight

DW = dry weight

1) Comment on the ATRO display mode

If the current measured value in the ATRO display mode is greater or less than the predefined limit value (i.e. greater than 999.99 %AD or less than –999.99 %AM), the ATRO result values are limited to 999.99%.

a – Weight in Grams


The weight of the sample is displayed (and printed out) in grams. With this setting, the Moisture Analyzer is used as a precision balance.

During the measurement current weight is constantly displayed in grams.

-%MC - Moisture Content

The moisture content of the sample is displayed (and printed out) as a percentage of the wet weight ($WW = initial\ weight = 100\%$).

During the measurement the value is constantly displayed in percent and as a graphical drying curve. The measured value is marked by "-%MC" (Moisture Content, e.g. -11.35 -%MC) also for the printed results and shown as a negative value.

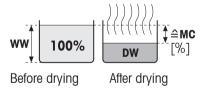

35

$$MC = -\frac{WW - DW}{WW} \cdot 100 \%$$

MC = Moisture Content [0...100%]

WW = wet weight

DW = dry weight

7.1.1.1.3 Start weight

Navigation: Home > Method Definition > Method name > Main measurement parameters > Start weight

The start weight affects both the duration of the measurement as well as the accuracy of the results. A small weight results in a shorter duration of the measurement but will decrease the accuracy of the result. With this function you can define a start weight so that the samples are approximately of equal weight thus improving the repeatability of the measurements. The weighing-in aid assists you in weighing-in the sample. At most of the samples the target weight is in the range of 2-5 g. We recommend to cover the whole surface of the sample pan with a thin uniform layer of sample.

The following parameters can be defined:

Parameter	Description	Values
Start weight	Defines a start weight so that the samples always have approximately the same weight.	OFF* I ON (0.5 g 101.090 g)
Weigh-in aid	Defines monitoring of target weight and tolerance (tolerance range: ±10%).	Passive* I Active
	Passive = tolerance is displayed. Active = tolerance is monitored. If the start weight is out of tolerance, the measurement cannot be started.	

^{*} Factory setting

7.1.1.2 Result and value handling

Navigation: Home > Method Definition > Method name > Result and value handling

The following parameters can be defined:

Menu item	Explanation	Further Information
Free factor	Defines a method-specific multiplication factor to calculate a corrected final result.	see [Free factor ▶ Page 36]

7.1.1.2.1 Free factor

Navigation: Home > Method Definition > Method name > Result and value handling > Free factor

This function allows you to multiply the result by a method-specific factor in the selected display mode. Furthermore, it allows you changing the result by an offset entry in the unit of the selected display mode. It is used to calculate a corrected final result (e.g., in order to compensate for systematic deviations from the reference result).



Note

- Free factor is not available in display mode "g" (weight).
- The calculation is carried out continuously during the measurement and displayed with the graphical drying curve.
- With free factor, the results are calculated and marked with an asterisk in the display.

The following parameters can be defined:

Parameter	Description	Values
Free factor	Defines the settings for free factor.	OFF* I ON
	OFF = no free factor applied.	
Factor	Defines the multiplication factor.	-10.000+10.000 (1.000)*
Offset	Defines the offset value in the selected display mode.	-1000.000 +1000.000 (0.000)*
Format	Defines the number of decimal places of the calculated final results, for displaying and printing.	x x.x x .xx x.xxx*

^{*} Factory setting

36

7.1.1.3 Workflow handling

Navigation: Home > Method Definition > Method name > Workflow handling

7.1.1.3.1 Starting mode

Navigation: Home > Method Definition > Method name > Workflow handling > Starting mode

In this menu you can choose whether the sample chamber performs automatic or manual functions (e.g. tare, stop drying).

The following parameters can be defined:

Parameter	Description	Values
Starting mode	Defines how the sample chamber is operated.	Automatic* I
		Manual

^{*} Factory setting

Automatic

Your instrument is set in the factory to the automatic operating mode. This mode can be used for most kind of samples. When closing the sample chamber, the weight of the sample is recorded and the measurement is started.

Manual

We advise you to use the manual operating mode for samples which contain readily volatile substances. The initial weight (wet weight) important for the determination of the moisture content is recorded. In the manual operating mode you have time for further preparation of the sample (e.g. mixing with quartz sand or even distribution of the sample) while weight losses due to evaporation during the preparation time are measured from the outset. As soon as the sample is ready for drying, close the sample chamber. As soon as the sample chamber is closed, drying starts. In the manual operating mode you can open the sample chamber during a drying operation. In contrast to the automatic mode, drying will not be stopped but simply interrupted until the sample chamber is closed again.

7.1.1.3.2 Accessories

Navigation: Home > Method Definition > Method name > Workflow handling > Accessories

For a method, certain accessories can be specified that must be used for the measurements. During the measurement, these accessories are then requested on the work screen.

The following parameters can be defined:

Parameter	Description	Values
Accessories	Defines the main accessory (pan) in which the sample is prepared.	Sample pan* I HA-Cage
Additional accessories (only available if the sample pan is the main accessory:	Defines additional accessories to be used with the sample pan.	☐ (deactivated)* I ☐ (activated)
 Glass fibre filter 		
 Second glass fibre filter 		
 Downholder 		

^{*} Factory setting

7.1.1.4 General method properties

Navigation: Home > Method Definition > Method name > General method properties

Moisture Analyzer Software Description

7.1.1.4.1 Method name

Navigation: Home > Method Definition > Method name > General method properties > Method name The following parameters can be defined:

Parameter	Description	Values
Method name	Renaming a method. Name should be unique and unambiguous.	any

7.1.1.4.2 Sample preparation

Navigation: Home > Method Definition > Method name > General method properties > Sample preparation

If a method is created with the method wizard (see [Creating a method with the Method Wizard ▶ Page 38]), the wizard automatically fills the **Sample preparation** field with the recommended preparation steps. The recommended preparation can be edited by tapping **Sample preparation**, but the number of characters in the text is limited to 100.

The sample preparation recommendation is not visible in the workflow during routine measurements.

The following parameters can be defined:

Parameter	Description	Values
	Specifies the sample preparation for a method recommended by the Method Wizard.	any

7.1.2 Creating a method

Prerequisite

The parameter menu appears only if a method already exists and is selected in the method definition dialog or if a new method has been created.

Main menu	Submenu	Further information
Main measurement parameters	Drying program (including temperature and switch-off criteria)	see Settings of Drying Program
	Display mode	see Settings of Display Mode
	Start weight	see Settings of Start Weight
Result and value handling	Control limits	see Control limits
	Resolution	see Resolution
	Free factor	see [Free factor ▶ Page 36]
	QuickPredict	see QuickPredict
Workflow handling	Starting mode	see Starting Mode
	Accessories	see [Accessories ▶ Page 37]
General method properties	Method name	see [Method name ▶ Page 38]
	Sample preparation	see [Sample preparation ▶ Page 38]

7.1.2.1 Creating a method with the Method Wizard

Introduction

The Method Wizard is a feature that offers assistance in developing a suitable moisture method for specific samples. Based on the characteristics of your sample as well as a test of the drying temperature and switch-off criterion, the Method Wizard is able to formulate a suitable method for your sample.


The Method Wizard guides you through the development of your moisture method in three simple steps:

- 1. **Questionnaire**: proposal of method parameters based on sample characteristics.
- 2. **Temperature assistant**: determination of drying temperature.

3. **Test measurement**: verification of temperature and selection of appropriate switch-off criterion After each step of the Method Wizard you can decide whether to accept the current method settings or to complete another step of the wizard. For steps 2 and 3, a reference moisture content value is needed.

Creating a new method with the Method Wizard

- Reference value (for Steps 2 and 3) is available.
- Method definition is selected.
- 1 Tap [**New...**].
- 2 Tap Define using Method Wizard.
 - The Method Wizard disclaimer text appears.
- 3 Read the disclaimer text and confirm by tapping [Next].
 - The keypad appears.
- 4 Enter a name for the new method. The name should be unique and unambiguous. 1 to 30 characters are possible (including spaces).
- 5 Confirm with [OK].
 - A short introduction appears.
- 6 Read the instruction and tap **Next**.
 - The Method Wizard starts.
- 7 Step 1 (Questionnaire): Follow through the questions by tapping the answers and then tapping [Next].
 - After all questions are answered, you are proposed a method and asked if you want to adopt this method or continue with the next step.
- To continue, tap Continue with temperature assistant (recommended).
 - To leave the wizard, tap Accept method parameters and exit wizard.
- 9 Confirm by tapping [Next].
- 10 **Step 2 (Temperature assistant):** Follow the instructions of the wizard to determine a suitable drying temperature. For questions about the temperature assistant, **see** Temperature assistant. This step may require several iterations.
 - After successful testing, you are proposed a drying temperature and asked if you want to complete the next step or adopt the current method.
- 11 To continue, tap Continue with test measurement (recommended).
 - To exit the wizard, tap Accept method parameters and exit wizard.
- 12 Tap [Next].
- 13 **Step 3 (Test measurement):** Follow the instructions of the wizard to select an appropriate switch-off criterion. For questions about the test measurement, **see** Test measurement.
 - After successful testing, you are recommended a switch-off criterion and are asked if you want to adopt it or perform another test.
- 14 To adopt the method, tap the desired switch-off criterion.
 - To reject the proposed switch-off criterion and perform another test at a different drying temperature, tap **Start over again using a different drying temperature**.
- 15 Confirm by tapping [Next].
- 16 To store the method, tap [Finish].
 - The method settings appear. The new parameters are already filled in.
- 17 **NOTICE: Data loss:** To definitively store your method settings, tap [**Save**] before you exit the method settings. Otherwise all data is lost.

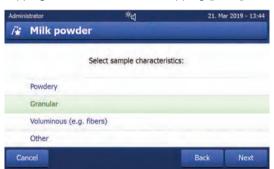

39

7.1.2.2 Creating a method manually

- Method definition is selected.
- 1 Tap [**New...**].
- 2 Tap Define method manually.
 - → The keypad appears.
- 3 Enter a name for the new method. The name should be unique (a method name can only exist once) and unambiguous. 1 to 30 characters are possible (including spaces).
- 4 Confirm with [OK].
 - → The parameter menu for the new method appears.
- 5 Set the desired parameters e.g. Main measurement parameters.
- 6 To store the method, tap [Save].

7.1.2.3 Creating a method by consulting the method library

The online METTLER TOLEDO moisture method library offers a wide range of tested and validated methods for different industries and substances. Optimize your moisture analysis with the help of our moisture expertise and profit from a wealth of know-how and widespread support.

Find the moisture library under:

www.mt.com/moisture-methods

Downloading a method from the method library

- PC with internet connection is available.
- Storage device (e.g. USB Stick) with at least 500 MB memory space is available.
- 1 Open the method library (see above) on your PC.
- 2 Click Moisture Applications.
- 3 Fill out the form according to the instructions to find a moisture method for your sample.
- 4 Click the method you wish to download.
 - → The access form opens.
- 5 Fill out the form and click **Download** to download the method file.
- 6 Transfer the file onto your storage device.

Importing the method to the istrument

- Method Definition is opened.
- 1 Connect your storage device to the instrument.
- 2 Tap [Import].
 - → Import methods screen opens.
- 3 Tap [Import from file].
 - → The list of connected hard drives opens.
- 4 Tap the device and location where your method file is stored and select your file.
- 5 Confirm by tapping **Import**.
- Your new method is imported and is now ready for routine measurements in the method list.

7.1.3 Editing a method

Editing an existing method

- Method definition is selected.
- 1 Tap the method in the list you want to edit.
 - → The parameter menu for the selected method appears.
- 2 Edit the desired parameters e.g. **Main measurement parameters**.

3 To store the method, tap [Save].

Copying an existing method

- Method definition is selected.
- 1 Tap the method in the list you want to copy.
 - The parameter menu of the desired method appears.
- 2 Tap [**Save as...**].
 - → The keypad appears.
- 3 Enter a new name for the copied method. The name should be unique and unambiguous. 1 to 30 characters are possible.
- 4 To store the method, tap [Save].

Removing an existing method

Note

All results of this method (incl. results of previous versions) as well as corresponding shortcuts are also removed.

- Method definition is selected.
- 1 Tap the method in the list you want to remove.
 - The parameter menu of the desired method appears.
- 2 Tap [Delete].
 - A message box appears.
- 3 Confirm with [Delete].
- → The desired method is removed.

7.2 Measurement

Navigation: Home > Measurement

Introduction

Before a measurement can be performed, a method first needs to be defined. To define a suitable method for your sample, **see** [Method definition > Page 30].

The instrument guides you through the steps during a measurement and displays results in real time. Nevertheless, please note the instructions in this manual, especially at the beginning of your work with the Moisture Analyzer, as a correct execution of the work steps and the handling of the samples can have a great influence on the results.

7.2.1 Performing a measurement

Before performing a measurement, a method has first to be defined. To find a suitable method for your sample, see Method Definition.

- The instrument must be connected to the power supply for approximately 60 minutes in order to reach operating conditions.
- To switch on the Instrument, press ♂.

Selecting a method

- 1 Tap on **Measurement**.
 - The methods list is displayed.
- 2 Select the method to determine your sample.
 - → The work screen of the method appears.
- 3 Open the sample chamber.


Moisture Analyzer Software Description

Placing the accessories

- The display now prompts you to load the predefined accessories and tare the balance.
- 1 Place the empty accessories in the sample pan handler.
- 2 Place the sample pan handler in the sample chamber. Ensure that the tongue of the pan handler fits exactly in the slot of the draft shield element. The sample pan must lie flat in the pan holder.

Note

We advise you to work with the sample pan handler at all times. The pan handler is ergonomic, automatic positioning, safe and provides protection against possible burns due to the hot accessories.

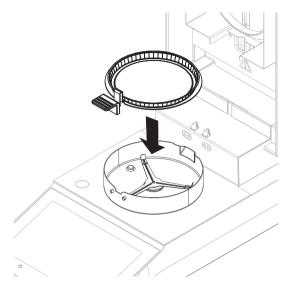

Taring the balance




Note

Before the sample pan is placed into the sample chamber, the left corner of the display must show **Open cover**, **place sample pan and tare**.

If the display shows Add sample, tap Cancel before placing the empty sample pan in the sample chamber.

- 1 Close the sample chamber.
 - → The instrument tares the balance (Starting mode: Automatic).
- 2 After tarring, open the sample chamber.
- After taring, the display prompts you to add the sample to the sample pan.
- 1 Add the sample to the sample pan. If you have defined a start weight, weigh the sample using the weighing-in aid.
- 2 Close the sample chamber.
- 3 Tap Start drying.
 - → The drying process starts.

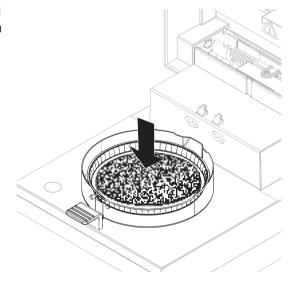

Drying process

The measurement process can be followed on the display.

- The drying process is continuously displayed graphically.
- The current temperature in the heating module is displayed as well as the elapsed drying time and the current drying value.
- The display shows the selected settings.
- The drying process can be aborted by tapping on Stop drying.

At the end of the drying process you can read the moisture content of your sample on the display.

The maximal drying time is 480 minutes (8 hours).

Removing sample




CAUTION

Burns due to hot surfaces!

Parts of the instrument can reach temperatures that cause injuries if touched. Sample, sample pan and other parts inside the sample chamber may still be hot.

- Do not touch the place marked with the warning symbol.
- The drying process is completed.
- Open the sample chamber.
- 1 Carefully remove the sample pan handler from the sample chamber.
- To perform a further measurement with the current method, tap Next Sample.
 - To perform a measurement with a new method, tap **Measurement**.
 - To return to the home screen, press 🔒.

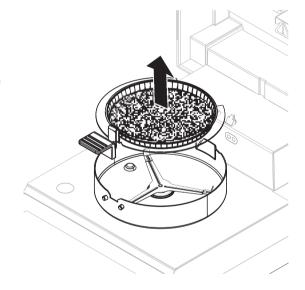

43

Stop drying

When you stop the measuring process by tapping [Stop drying], you can choose between two options:

- Abort measurement and save data
 - Already sampled data are stored and an entry will be made in the results. The result is marked as aborted.
- Cancel

The process will be continued.

Adding a comment

At the end of a measurement you can add a comment to the result of the measurement. This comment is stored in the measurement result and can be printed. A comment can only be entered before leaving the current measurement.

- 1 To create a comment, tap **Note**.
 - The keyboard dialog appears.

- 2 Enter a comment.
- 3 Confirm with **OK**.

7.2.2 Optimum sample preparation

Preparation of the sample is decisive for the speed of the measurement process and the quality of the measurement results.


Basic rules for the preparation of samples:

- The amount of sample you select should be as small as possible but as large as necessary.
- The more inhomogeneous the sample substance, the larger the amount of sample needed to obtain a repeatable result.
- Distribute the sample evenly over the sample pan. This increases the surface area of the sample and facilitates heat absorption.
- With liquid, fat-containing, melting, highly reflecting samples or samples which form a skin on their surface
 when heated, use the sample with the glass fiber filter available as optional equipment, see Accessories
 and Spare Parts.

7.2.3 Work with weighing-in aid

The weighing-in aid can be defined for each method and facilitates weighing in of the sample to a target value. This is particularly necessary if you require all the samples of a method you wish to process to have same weight in order to improve the repeatability of the measurement results. In addition, the weighing-in aid can be set as active so that the drying process cannot be started if the weight of sample is outside the set tolerance. You are therefore compelled to weigh-in the correct quantity of sample. If all sample weighed are within tolerances then this will improve the repeatability. The weighing-in aid is available only if start weight has been activated.

For further information **see** [Settings of Start Weight ▶ Page 36].

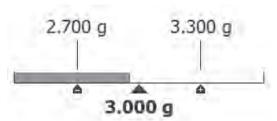

Icon	Function
	Lower weight limit (tolerance range)
	Target weight
△	Upper weight limit (tolerance range)

7.2.4 Working with shortcuts

Navigation: Home > Measurement > Method name > 🗗

With shortcuts, methods can directly be started from the home screen. Shortcuts are being created for important methods that are often used.

Create a shortcut

- 1 Tap on **Measurement**.
 - The list of the defined methods appears.
- 2 Tap on **Edit shortcuts...** that can be found in the action bar.
- 3 Tap on the method from the list for which you want to create the shortcut.
 - → The window **Shortcut name** appears.

4 Enter a shortcut name.

Note

The shortcut name has a maximum length of 30 characters. Two lines with about 16-20 characters of the shortcut name are being shown in the shortcut on the home screen. The space button can be used to break the line. We recommend to use not more than 10 characters per line.

- 5 Tap on **OK**.
- 6 Tap on Save.
- → The new shortcut appears on the home screen.

Editing a shortcut

- 1 Tap on **Measurement**.
 - → The list of the defined methods appears.
- 2 Tap on Edit shortcuts....
- 3 Tap on the defined method from the list for which you want to edit the shortcut.
- 4 Define the shortcut name.
- 5 Tap on **OK**.
- 6 Tap on Save.
- The shortcut has been edited.

Removing a shortcut

- 1 Tap on **Measurement**.
 - The list of the defined methods appears.
- 2 Tap on Edit shortcuts....
- 3 Disable the shortcut that should be removed from the home screen.
- 4 Tap on Save.
- → The shortcut is removed from the home screen.

7.3 Results

Navigation: Home > Results

The device can save a total of 100 measurement results. When there are more than 100 results, the oldest result will automatically be overwritten. Therefore we recommend to safe the results frequently.


7.3.1 Graphical evaluations of measurement results

Navigation: Home > Results

Moisture Analyzer Software Description

This function allows you to manage and evaluate measurement results.

Method list

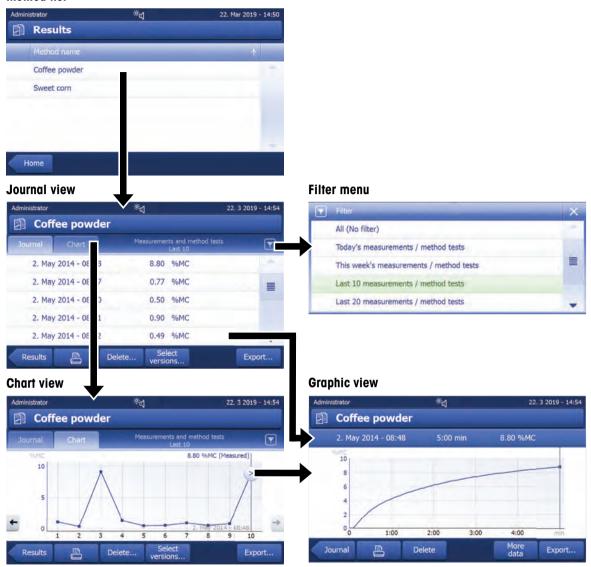

1 Method list

- Tap the method you want to evaluate.
 - The journal view appears.

2 Journal view

46

The journal view allows you to start different graphical evaluations from a measurement series. You can perform the following functions:

- To start filter function, tap \(\bigvee\).
 - The filter menu appears.
- To start chart view, tap Chart.
 - → The chart view of the series of measurement appears (4).
- To start graphic view, tap a measurement result.
 - The graphic view appears (5).

3 Filter menu

The filter function allows you to evaluate the measurement series according to various criteria. You can select the following criteria:

- All (No filter)
- Today's measurements
- This week's measurements
- Last 10 measurements*
- Last 20 measurements
- Advanced filters
- * Factory setting
- Tap **X** to close filter menu.

Advanced filters

The menu **Advanced filters** offers you a number of additional filter criteria.

The following settings can be defined:

Parameter	Description	Values
Time range	With the option Date range the results are being filtered by date. Begin and end date must be defined.	No time range limit I Date range I Last
	With the option Last measurements the last measurements and method tests that have been performed can be shown. The	measurements
	number must be between 1 and 100.	

4 Chart view

This view allows you to display the results of a series of measurements depending on the filter settings. The narked measuring point is displayed with date, time, and the measurement result. You can perform following functions:

- To jump forward to the next measurement result, tap -> .
- To jump backwards to the previous measurement result, tap <- . Alternatively, tap directly on the desired
 measurement point.
- To recall the full measurement curve of the corresponding measurement, tap 2.

Moisture Analyzer Software Description

5 Graphics view

With this function you can display a graphical view of the detailed results of a single measurement.

You can perform the following functions:

Printing result

To print the result, tap \(\bullet \).

Exporting result

To export the result, tap Export.

Deleting result

- To delete this measurement result, tap **Delete**.

With the auto-delete functionality the oldest results will automatically be removed.

More data

To display more measurement data, tap More data

With this function you can display a graphical view of the detailed results of a single measurement.

7.3.2 Method versioning

Navigation: Home > Results

Every method has a method versioning function that works in the background. Whenever the settings of a method are changed, the versioning function automatically generates a new version of the method. This has also an effect on the results: Every measurement overtakes the current version of the method.

Measurement and method test results can be filtered based on the method version as follows:

- 1 Open the menu Results.
- 2 Select a method.
 - → A journal with the results of the performed measurements and method tests appears. If there are results in several versions, the button Select versions... appears in the action bar.
- 3 Tap Select versions....
 - → The dialogue Method versions appears.
- 4 Select the versions you want to filter.
- 5 Tap **Save** to confirm the selected versions.
- The system returns to the result list. The results based on the selected versions appear on the list, the other results are being grayed out.



The results filtered by the method version can be exported or printed. If the results are being exported as PDF or being printed, the results of the different versions are being listed based on their version.

7.3.3 Exporting results

7.3.3.1 Exporting results to a storage device

Navigation: Home > Results

Results can be exported to an external storage device e.g. a memory stick. Data in CSV format can be imported into e.g. MS Excel for further evaluations.



Note

Results cannot be imported in a Moisture Analyzer.

Sample procedure

- Result is activated.
- An external storage device is connected e.g. memory stick.
- 1 To start, tap **Export**

- **Export results** appears.
- 2 Tap Export target > Peripheral
- 3 Tap Location > Browse and select.
 - Location appears.
- 4 Select location for file and confirm with **OK**.
- 5 Enter new file name, if needed.
- 6 To start export, tap Export.

7.3.3.2 Exporting results with EasyDirect Moisture

With **EasyDirect Moisture**, measurement and routine testing results are collected in a database. Results can be viewed, exported and also routine tests can be scheduled. The PC software must be installed on an external PC and the device must be connected to that PC with a USB cable or RS232 cable.

Before the function **EasyDirect Moisture** can be used, the following settings must be defined on the instrument:

Defining EasyDirect Moisture as peripheral device

Navigation: Settings > Instrument settings > Peripherals.

- Peripherals is opened.
- No printer is installed as a peripheral. Uninstall printer if necessary.
- 1 Tap [**New...**].
 - → The list of possible peripherals appears.
- 2 Select [EasyDirect Moisture].
 - → The parameters for EasyDirect Moisture appear.
- 3 Under **Device interface**, select the desired interface for the EasyDirect Moisture connection:
 - For a RS232-interface, select RS232.
 - For a USB-interface, select USB.
- 4 Set the rest of the settings.
- 5 To save the settings, tap [Save].
- → The option EasyDirect Moisture has been defined as peripheral device on the instrument.

See also

Peripherals ▶ Page 61

7.3.3.3 Export settings

7.3.3.3.1 Single result export settings

The following parameters can be defined:

Parameter	Description	Values
File name	Defines the name of the new data file.	any
File type	Defines the type of the data file.	CSV* I PDF
Language	Defines the language for export files.	English I Deutsch I Italiano I Français I Español I Português I Polski I Magyar I Český I Русский I 日本語 I 中文 I 한국어/조선말
Export target	Defines the destination of data transfer.	Peripheral

^{*} Factory setting

Moisture Analyzer Software Description

Peripheral

Parameter	Description	Values
Location	Finds the location on the storage device.	Browse and select

7.3.3.3.2 Multiple result export settings

Summary export

The summary export contains method parameters and the end results per measurement. The intermediate values are not exported.

Multiple export

- The multiple export contains method parameters, method data, intermediate values and end results per measurement.
- Each measurement is exported as a single file.

The following parameters can be defined:

Parameter	Description	Values
File name prefix	Defines the name of the new data file.	any
Export selection	Custom selection of multiple results for the export from selected measurements. The selection can be made via the filter function.	Selection: ☐ I ✓
Export content	Defines the content of the multiple results export.	Summary export I
	Summary export (one file, no intermediate values): Results as summary report in one file. Multiple export (one file per measurement): Results with intermediate measurement values as separate files with the same layout as for the single result export.	Multiple export
File type	Defines the type of the data file.	CSV* I PDF
Language	Defines the language for export files.	English I Deutsch I Italiano I Français I Español I Português I Polski I Magyar I Český I Русский I 日本語 I 中文 I 한국어/조선말
Export target	Defines the destination of data transfer.	Peripheral

Peripheral

Parameter	Description	Values
Location	Finds the location on the storage device.	Browse and select

^{*} Factory setting

7.4 Test/Adjust

50

Navigation: Home > Test/Adjust

Introduction

For optimum precision and repeatable results of the Moisture Analyzer, the function of the instrument should be tested on a regular basis and, if necessary, adjusted. The frequency of testing and adjusting depends on the use and location of the instrument.

Main menu	Submenu	Further information
Adjustments	Weight adjustment - internal see Weight Adjustment - Internal	
	Weight adjustment - external	see Weight Adjustment - External
	Temperature adjustment see Temperature Adju	

Tests	Weight test - internal	see Weight test - internal
	Weight test - external	see Weight test - external
	Weight test - Repeatability external	see Weight test - Repeatability external
	Temperature test	see Temperature test
	SmartCal x1 test	see [SmartCal test ▶ Page 53]
History	Weight test history	see History
	Weight adjustment history	
	Temperature adjustment history	
	Temperature test history	
	SmartCal x1 test history	

7.4.1 Tests

Navigation: Home > Test/Adjust > Tests

In this menu item you can check several settings, such as the adjustment of the balance or the temperature of the heating module. For relevant settings **see** [Test / adjustment settings \blacktriangleright Page 66].

Introduction

Routine tests are performed to ensure accurate results. Therefore the tests should be performed in fixed, regular intervals depending on your Quality Management system and the results should be documented in a traceable way.

Based on your process requirements, METTLER TOLEDO can help you to define the routine tests to be performed. For further information, contact your local METTLER TOLEDO representative.

7.4.1.1 Temperature test

Navigation: Home > Test/Adjust > Tests > Temperature test

Target of the test

The temperature test checks if the temperature sensor is working correctly and if it delivers correct results. Using the temperature kit, the difference of the measured temperature at the kit (target temperature 1) and the sensor (target temperature 2) is used to calculate the precision of the sensor.

It is possible to start a test and subsequently transform it into an adjustment, if required.

Equipment

• Temperature kit.



CAUTION

Burns due to hot surfaces!


Parts of the instrument can reach temperatures that cause injuries if touched.

- 1 Let the instrument cool down before setting up.
- 2 Hold the temperature kit by its handle only.


Moisture Analyzer Software Description

Installing the temperature kit

- Instrument is cooled down to < 50 °C.</p>
- Instrument is switched on.
- Sample chamber is open.
- 1 Remove sample pan handler (with pan) (1).
- 2 Remove sample pan holder (2).
- 3 Clean the contact areas (3) if needed.

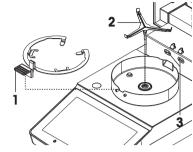

4 Insert Temperature kit (4).

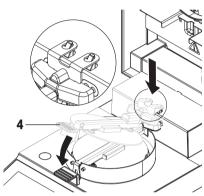

Testing procedure

- Test settings are defined, **see** Test Settings.
- Temperature kit is installed.
- 1 Tap Temperature test.
 - → Work screen Temperature test appears.
- 2 Close the sample chamber.
- 3 Tap [Start test].
 - → The testing process starts: The instrument first heats up to target temperature 1 (lower) for 15 minutes, then, for another 15 minutes, heats further up to target temperature 2 (higher). The screen shows the target temperature values (Target), measured at the sensor, and the actual temperature (Actual), measured at the temperature kit.
 - → The instrument indicates whether the test has passed or failed according to defined tolerances.
- 4 If an adjustment is needed, decide if you want to perform it now, see Temperature adjustment.
- To return to the menu Tests, tap [Tests].
 - To print the results, tap [].
 - To display or print the results at any time, tap [History].
 - To return to the home screen, press [1].
- 6 Let the instrument cool down before removing the temperature kit.

See also

52

- Accessories and Spare Parts ▶ Page 87
- Test settings ▶ Page 67
- Temperature adjustment ▶ Page 56
- Accessories ▶ Page 87

7.4.1.2 Weight test - external

Navigation: Home > Test/Adjust > Tests > Weight test - external

Target of the test

The external weight test determines if the integrated balance and the weight value display are working correctly by placing a test weight. The test ensures correct and precise results.


Equipment

Test weights. To guarantee full traceability, METTLER TOLEDO recommends to use certified weights.

Testing procedure

- Test weight and test settings are defined.
- 1 Open sample chamber.
- 2 Tap the item Weight test external.
 - → The test screen appears.
- 3 Clear sample pan holder and tap Start test.
 - The required value flashes and prompts you to load the weight.
- 4 Load the required test weight.
 - The balance indicates the test process, ——— appears.
- 5 When the display flashes **000.00** remove the test weight.
- The display shows the test results and indicates whether the test has passed or failed according to defined tolerances.

To print the results, tap [具].

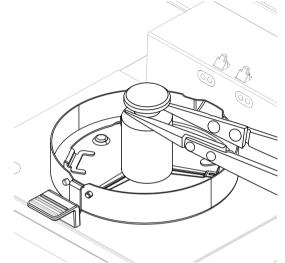

53

Timeout

- In case of a timeout of over 75 seconds, the error message appears Weight out of range. The test has been aborted. The test was not executed.
- 1 Confirm with [OK].
- 2 Repeat the test.

See also

Accessories ▶ Page 87

7.4.1.3 SmartCal test

Target of the test

A SmartCal test is a quick and integral test of the instrument functions. SmartCalTM is a granular test substance whose moisture content value, when heated at a specific temperature for a specific time, is precisely defined. By performing a measurement with SmartCalTM and comparing the results to the control limits, it can be tested if both the integrated balance as well as the temperature sensor work well together and if the instrument functions correctly.

However, SmartCal tests do not supersede the recommended periodic weight or temperature tests.

There are four SmartCal tests with four pre-defined temperatures and two additional tests where temperature and control limits can be defined by the customer.

Equipment

- SmartCal[™] Moisture Analyzer Test Substance.
- For best results and easy handling, use a RHT sensor.
 If no RHT sensor is available, the room temperature and relative humidity must be entered manually.

Certified testo RHT sensors are supported.

Certified RHT sensors

The following RHT models from Testo are supported: testo 435-2, testo 435-4, testo 635-2, testo 735-2.

The testo RHT sensor has to be connected to the USB port on the terminal.

For more information about Testo, see

http://www.testo.com

Rules for handling SmartCal™

- Store SmartCal[™] at room temperature.
- Remove SmartCal[™]-stick from the blister pack right before use.
- Do not use damaged or expired sticks (the expiry date is printed on the blister pack and on the stick (e.g. Exp08.2021).
- After a SmartCal[™] test, the test substance can be disposed of as normal waste.

Additional information about SmartCal™ can be found under:

www.mt.com/smartcal

Testing procedure

The test is carried out in the same way as a measurement.

Note

Every SmartCal test has a maximum number of 30 test results. If there are more than 30 results for one test, the oldest results will automatically be deleted.

- The instrument has been connected to the power supply for at least 60 minutes in order to reach operating conditions.
- If available, a RHT sensor is connected.
- The SmartCal package is acclimatized.
- Test temperature and control limits are set.
- The sample chamber is open.
- 1 Tap [SmartCal x1 test].
 - The method list is shown on display.
- 2 Select a method from the list.
 - → The work screen appears.
- 3 If the test is performed without RHT sensor, tap [Enter room temperature and humidity], enter current room temperature and humidity and confirm with [OK].
- 4 Enter SmartCal[™] lot number if needed. Tap [**OK**]
- 5 Place the sample pan handler with the aluminum sample pan in the instrument.
- 6 Tap [->0/T<-] to tare the balance.
- 7 Remove a SmartCal[™] stick from the blister pack, open it and distribute the entire contents evenly over the sample pan. If necessary, carefully rotate and tilt the sample pan until it is fully and uniformly covered with granulate.
- 8 Start the test immediately by tapping [Start drying].
 - The SmartCal test procedure starts.
 - ➤ When the test is finished, the instrument indicates whether the test has passed or failed. The display shows the normalized result. If the SmartCal result has failed, see [Measures after failed test ▶ Page 55].
- To return to the menu Tests, tap [Tests].
 - To print the results, tap [].
 - To display or print the results at any time, tap [History].
 - To return to the home screen, press [1].

See also

- Test settings ▶ Page 67
- Accessories ▶ Page 87

7.4.1.4 Measures after failed test

In case of a failed test, take the following measures:

- 1 Repeat the test and ensure that all steps have been carried out correctly.
- 2 If the failure persists, perform an according adjustment.
 - If a weight test has failed, perform a weight adjustment.
 - If a temperature test has failed, perform a temperature adjustment.
 - If a SmartCal test has failed, perform a weight adjustment first, then perform another SmartCal test and, if the failure persists, perform a temperature adjustment.
- 3 If the failure persists, **see** Malfunctions when operating the instrument.
- 4 If the failure persists, contact your METTLER TOLEDO service representative.

7.4.2 Adjustments

Navigation: Home > Test/Adjust > Adjustments

Introduction

To guarantee correct results and high performance of the instrument, it is necessary to adjust its functions:

- After setting up the balance for the first time.
- After a change of location.
- If a test indicates that an adjustment is due.

Test the instrument exclusively under operating conditions.

METTLER TOLEDO offers an adjustment service – please contact your METTLER TOLEDO representative.

See also

Test / adjustment settings ▶ Page 66

7.4.2.1 Weight adjustment - external

Navigation: Home > Test/Adjust > Adjustments > Weight adjustment - external

Target of the adjustment

In an external weight adjustment, the integrated balance is adjusted by placing a predefined weight as a reference for the instrument to re-calibrate itself.

Equipment

Test weights. To guarantee full traceability, METTLER TOLEDO recommends to use certified weights.


Testing procedure

- Test weight and test settings are defined.
- 1 Open sample chamber.

Moisture Analyzer Software Description

- 2 Tap the item Weight test external.
 - → The test screen appears.
- 3 Clear sample pan holder and tap Start test.
 - The required value flashes and prompts you to load the weight.
- 4 Load the required test weight.
 - → The balance indicates the test process, ———— appears.
- 5 When the display flashes **000.00** remove the test weight.
- The display shows the test results and indicates whether the test has passed or failed according to defined tolerances.

To print the results, tap [].

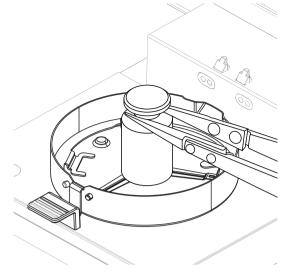

Timeout

- In case of a timeout of over 75 seconds, the error message appears Weight out of range. Adjustment aborted. The test was not executed.
- 1 Confirm with [OK].
- 2 Repeat the adjustment.

See also

Accessories ▶ Page 87

7.4.2.2 Temperature adjustment

Navigation: Home > Test/Adjust > Adjustments > Temperature adjustment

This function allows you to adjust the temperature control of the heating module. You need the optional Temperature Kit for performing this function, **see** [Accessories and Spare Parts ▶ Page 87]. To learn when an adjustment of the heating module is necessary, **see** Notes on adjustment of balance and heating module. We advise you to wait at least 30 minutes after a drying operation (or previous heating module adjustment) before performing the adjustment.



56

Note

- The duration of the temperature adjustment is 30 minutes. (15 minutes for each measuring temperature).
- The sample chamber must be less than 50 °C before a further adjustment or test can be done.
- It is also possible to perform a test first, and then to transform the test into an adjustment if required. Only
 possible for tests conducted with 2 temperatures, see [Temperature test ▶ Page 51].
- For temperature adjustment settings see [Adjustment settings ▶ Page 69].



CAUTION


Burns due to hot surfaces!

Parts of the instrument can reach temperatures that cause injuries if touched.


- 1 Let the instrument cool down before setting up.
- 2 Hold the temperature kit by its handle only.
- Adjustment settings are defined.
- Tap menu item Temperature adjustment.
 - → Work screen Temperature adjustment appears.

Installing the temperature kit

- Instrument is cooled down to < 50 °C.
- Instrument is switched on.
- Sample chamber is open.
- 1 Remove sample pan handler (with pan) (1).
- 2 Remove sample pan holder (2).
- 3 Clean the contact areas (3) if needed.

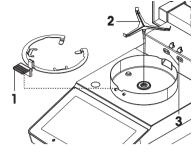

4 Insert Temperature kit (4).

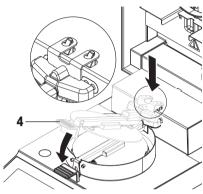

57

Start adjustment

- Draft shield element is installed.
- Sample pan handler and sample pan holder is removed.
- **Temperature kit** is inserted.
- 1 Close sample chamber.
- 2 Tap [Start adjustment].

The adjustment process takes 30 minutes.

- → The instrument starts the adjustment process. The screen shows the current temperature and the remaining time.
- → The instrument heats up to the target temperature 1 for determining the predefined lower temperature.
- After 15 minutes, the instrument adjusts the lower temperature and continues with the process.
- → The instrument heats up to the target temperature 2 for determining the predefined upper temperature.
- → After 15 minutes, the instrument adjusts the upper temperature and concludes the adjustment.
- To print out the results, tap \(\bullet
 - To return to the home screen, tap .
- 4 After cooling down, remove **Temperature kit** and set up drying unit for measurement.

See [History ▶ Page 58] and [Setting up the instrument ▶ Page 14].



Note

Your calibrated temperature kit can be re-calibrated. Please contact your local METTLER TOLEDO representative for more information.

7.4.3 History

Navigation: Home > Test/Adjust > History

The instrument always records all test or adjustment operations that have been made and saves them in a memory which is specially protected against power failure (only the last adjustment or test is being displayed).



Aborted adjustments or tests will not be saved.

When the History is selected, a list appears with the history selection as follows:

Weight test history

Temperature adjustment history

Weight adjustment history

SmartCal x1 test history

• Temperature test history

The following information is displayed:

Weight test history

Weight test external	
Calibration weight identification	
Nominal weight	g
Actual weight	g
Difference	g
Tolerance	g
Temperature	°C

Weight adjustments history

Weight adjustment external	
Calibration weight name	
Calibration weight identification	
Nominal weight	g
Temperature	°C

Temperature adjustment/test history

Temperature adjustment		Temperature test	
Adjustment kit identification		Adjustment kit identification	
Temperature 1	target - actual	Temperature 1	target - actual
Temperature 2	target - actual	Temperature 2	target - actual
		Free temperature	target - actual
		Tolerances for each temperature	°C

SmartCal x1 test history

Following parameters are available for each SmartCal method.

SmartCal test	
User name	
SmartCal identification	
Start weight	g
Total test time	minutes
Dry weight	%MCN
Test result	passed / failed
Date and time	
Leveling information	

The selected individual results can be printed out by tapping [].

7.5 Settings

Navigation: Home > Settings

In this section you will find information how to define the general settings of your instrument.

Main menu	Submenu	Further information	
User preferences	Languages	see [Languages ▶ Page 59]	
	Screen	see [Screen ▶ Page 60]	
	Sound	see [Sound ▶ Page 60]	
	Protection	see [Protection ▶ Page 60]	
Instrument settings	Regional Settings	see [Regional settings ▶ Page 61]	
	Peripherals	see [Peripherals ▶ Page 61]	
	Touch screen adjustment	see [Touch screen adjustment ▶ Page 63]	
Application settings	Identifications management	see Identifications management	
	Print and export management	see [Print and export management ▶ Page 64]	
Quality management	Test / Adjustment settings	see [Test / adjustment settings ▶ Page 66]	
System and data management	Export / Import	see Export / import	
	Back up / Restore	see [Back up / Restore ▶ Page 71]	
	Export event logs	see [Export event logs ▶ Page 72]	
	Reset	see Reset	
	Update	see Update	
Help and tutorials	Instrument tutorial	see Instrument tutorial	

7.5.1 User preferences

Navigation: Home > Settings > User preferences

7.5.1.1 Languages

Navigation: Home > Settings > User preferences > Languages

Moisture Analyzer Software Description

The following languages can be chosen:

Parameter	Description	Values
Display language	Defines user-specific display language.	English Deutsch Italiano Français Español Português Polski Magyar Český Русский 日本語 中文 한국 어/조선말
Keyboard language	Defines user-specific layout of the keypad.	English I Deutsch I Italiano I Français I Español I Português I Polski I Magyar I Český

7.5.1.2 Screen

Navigation: Home > Settings > User preferences > Screen

The following parameters can be defined:

Parameter	Description	Values
Brightness: x1 %	Defines user-specific brightness of the display.	10100 % (80 %)*
Color scheme	Defines user-specific display color scheme.	Blue* Green Orange Grey

^{*} Factory setting

7.5.1.3 Sound

Navigation: Home > Settings > User preferences > Sound

The following parameters can be defined:

Parameter	Description	Values
Sound	Activate / deactivate of audible feedback and adjust the Volume: x1 %.	OFF* I ON I
	Audible feedback: • when tapping a button.	0100 % (50 %)*
	upon measurement completed.	
	when an error message appears.	

^{*} Factory setting

7.5.1.4 Protection

Navigation: Home > Settings > User preferences > Protection

The following parameters can be defined:

Parameter	Description	Values
	With this function activated the settings are protected from	ON I OFF
settings	accidentally changes through the operator.	



Note

With the function Protection of settings activated, the following functions stay activated:

- Measurement (edit shortcuts and deleting shortcuts is inactive)
- Results (Delete results is inactive)

Tests

7.5.2 Instrument settings

Navigation: Home > Settings > Instrument settings

7.5.2.1 Regional settings

Navigation: Home > Settings > Instrument settings > Regional Settings

The following parameters can be defined:

Parameter	Description	Values
Date	Sets the date of the clock of the instrument.	Day I Month I Year
Time	Sets the clock time of the instrument.	Hours I Minutes
Date format	Defines the date format. D = Day M = Month Y = Year	D. MMM YYYY* I MMM D YYYY I DD.MM.YYYY I MM/DD/YYYY I YYYY/MM/DD I YYYY-MM-DD
Time format	Defines the time format. M = Minutes	24:MM* 12:MM 24.MM 12.MM
Daylight saving time	Activates or deactivates the daylight saving time.	OFF* I ON

^{*} Factory setting

Daylight saving time

Define the following parameters:

Parameter	Description	Values
Offset	Defines the offset relevant for the daylight saving time.	30120 min (60 min)*
Start	Defines the point in time from which daylight saving time is active.	Time I Day of the week I Week of the month I Month
End	Defines the point in time from which daylight saving time is deactivated.	Time I Day of the week I Week of the month I Month



Note

The date and time of the moisture analyzer is relevant. The date and time of a connected printer will be ignored.

7.5.2.2 Peripherals

Navigation: Home > Settings > Instrument settings > Peripherals

In this menu item you can configure one printer and one host connection.

The **MT-SICS** host connection (remote client) is enabled by default and only usable with the METTLER TOLEDO specific protocol MT-SICS.

MT-SICS commands and functions can be transmitted via the RS232 interface (COM1 port) or the USB device interface. The use of the USB device interface requires the installation of an USB device driver on the PC, **see** [Installing USB device driver > Page 17].

Commands and further information please refer to the Reference Manual MT-SICS Interface Commands for Halogen Moisture Analyzers downloadable from the internet.

►www.mt.com/moisture

Installing Printers / Computers

No printer is installed by default. To install a printer:

- The printer is connected.
- The printer is powered on.
- 1 Tap **New...**.
 - Peripheral type list appears.
- 2 Select device.

Note

Currently supported devices from METTLER TOLEDO with preset parameters:

- P-2x Strip Printer
- P-4x Strip Printer
- P-5x Strip Printer
- EasyDirect Moisture
- 3 Change the parameters only if you have to. The parameters are predefined and any changes need advanced knowledge.
- 4 To store the settings, tap **Save**.
- 5 To activate the interface, tap **ON**.
 - ⇒ Status changes from **Disabled** to **Enabled**.



- A printer or host connection can also be deleted.
- For strip printer settings see Recommended Printer Settings (Strip printer).
- For defining the printouts see [Print and export management ▶ Page 64].
- Only one printer can be installed. To install another printer, you have to delete the old printer first.
- If EasyDirect Moisture is connected, the icon \(\bullet\) can be used to perform a manual transfer.

Parameters

The following parameters can be defined:

Parameter	Description	Values
Device Name	Defines a name for the device.	any
Device interface	Defines the device interface.	RS232* I USB
Device Port	Defines the device port. COM1: RS232 / SOU: USB Device (Driver required) Note It is possible to use an USB-to-Serial adapter. If such an adapter is connected, these devices are shown as e.g.SOU1, SOU2 (SOU stands for serial over USB)	COM1* I SOU
Baudrate	Defines the speed of the data transmission (data transfer rate / Baud rate).	1200 2400 4800 9600* 19200 38400 57600 115200
DataBits	Defines the number of data bits.	7 8*
Parity	Defines the variant of parity bit.	None* Odd Even
StopBits	Defines the stop bits of the transmitted data.	1 StopBit* I 2 StopBits
Flow Control	Defines the type of flow control (Handshake).	None I Xon/Xoff*

^{*} Factory setting

62

USB devices

Connecting USB strip printers

Strip printers connected over USB or with USB adapters need the following settings:

Device interface: RS232CDevice Port: SOUx

• **Baudrate**: 19200



- A bar code reader can be connected to the USB port. With a connected bar code reader, bar codes can be added in a keyboard dialog.
- If the printer is not powered on when adding it to the peripherals, the device port SOUx might not be visible.

7.5.2.3 Touch screen adjustment

Navigation: Home > Settings > Instrument settings > Touch screen adjustment


If the alignment of the touch-sensitive areas of the screen correspond not exactly with the position of the buttons e.g. after a software update was performed, then you can solve this error by executing this function.



NOTICE

Damage to the touch screen due to pointed or sharp objects

- Operate the touch screen with your fingers.
- Menu Instrument settings is opened.
- 1 Choose Touch screen adjustment and tap Start.
 - → Adjustment screen with a target circle appears.
- 2 Tap the center of the target until it moves to the next position on the screen. Tap the center of the target as accurately as possible and make sure not to touch any other locations on the screen and the fix key areas.
- 3 Repeat this procedure until a message appears (New calibration settings have been measured).
- 4 Tap the screen to confirm or wait for 30 seconds for canceling (timeout).
 - A message box appears.
- 5 To store the setting, tap **OK**.

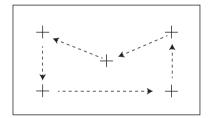

63

7.5.3 Applications settings

Navigation: Home > Settings > Application settings

7.5.3.1 Identifications management

Navigation: Home > Settings > Application settings > Identifications management

Identifications (ID) contain descriptive text for individual measurements, which enables samples to be easily allocated to specific tasks or customers. This feature allows you to define identifications in order to comment measurements, such as company id, id badge, id sample. In the ID panel on the work screen, the values (comments) can be entered or changed before measurement, **see** [Work screen \times Page 23]. IDs have an impact on all measurements, regardless of methods. The IDs are included on printouts or can be exported to a PC.

Main menu	Submenu	Further information
Identifications usage	Identifications input	see below
	Clear input after measurement	
Identifications (x1)	no submenu	

The following parameters can be defined:

Parameter	Description	Values
Identifications	Defines the method for entering comments.	OFF* I ON I
input	ON = activates ID panel on work screen.	Free input I
	Free input = the IDs appear after tapping on the ID panel (work screen) for entries or changes of values (comments).	Guided input
	Guided input = each ID is displayed individually in order of definition (ID1 to ID10) before measurement, to make entries or changes of values (comments).	
Clear input after measurement	Defines the ID whose value (comment) is deleted after measurement.	OFF* I ON I ID selection: ☐ I ☑ I
Identifications (x1)	Defines general identifications in order to comment measurements.	-l +*lany
	+ = New identifications (max. 10 IDs, 120 characters) - = Remove identifications (selectable)	

^{*} Factory setting

7.5.3.2 Print and export management

7.5.3.2.1 Print management

Print management (Strip)

Navigation: Home > Settings > Application settings > Print and export management > Print management (Strip)

If a strip printer is connected and activated, results can be automatically printed out. The intermediate values are printed out in the pre-selected intervals and the final results are logged at the completion of the measurement. The printout contains information to the instrument, time, date as well as the method and results, depending of the protocol. The level of detail depends on the selected type of printout (short, standard) and the activated protocol items (settings or method).



Note

- The definitions in this menu item are the default settings for the printouts. These parameters are displayed at every initialization of a Printout and can be modified again.
- . The setting for printout language in this menu item has no impact on test or adjustment printouts.
- Not all languages are compatible with some printers, and also depending on the settings of the printer, **see** Recommended printer settings (Strip printer).
- For installing and defining a printer see [Peripherals ▶ Page 61].

The following parameters can be defined:

Parameter	Description	Values
Printout type	Selects a detailed printout or in a short form.	Standard* Short
		Custom 1 Custom 2

Printout language	Defines the language for general printouts. Note No impact on test or adjustment printouts.	English I Deutsch I Italiano I Français I Español I Português I Polski I Magyar I Český
Print results automatically after measurement	Print results automatically after measurement if activated.	☐ (deactivated)* I ☑ (activated)
Include inter- mediate results to printout	Defines the time intervals at which intermediate results are transferred to the printer.	OFF* 10 s 30 s 1 min 2 min 10 min

^{*} Factory setting

Customized printouts

With the options **Custom 1** and **Custom 2** the information that appear on the printout can be selected individually. The information that appear on the printout can be defined in the sub-option **Printout content**. The following information for **Printout content** can be enabled/disabled or defined:

Parameter	Description	Values
Head	METTLER TOLEDO I Halogen Moisture Analyzer I Blank line	
S/N instrument and software	Type I SNR (Drying unit) I SNR (Terminal) I SW(Drying Unit) I SW(Terminal) I Blank line	
Adjustment info	Weight Adjustment I Temperature Adjustment I Blank line	
Method parameters	Method Name Drying Prog Drying Temp Switch Off Display m Weight Tol Weigh-in aid Start Mode Free factor Blank line	ode Start Weight Start
ID's	ID's I Blank line	
Measurement	Start weight Intermediate results Total time Dry Weight Moistu Offset Measured Result Measurement status Note Blank line	ıre Content Factor
Signature 1	☐ (deactivated)* I ☑ (activated)	
Signature 2	☐ (deactivated)* I ☑ (activated)	



Note

The information **Date / Time** and the end result are always activated and appear at the end of the printout. The settings for the option **Intermediate results** are still set but do only appear on short and standard printout.

7.5.3.2.2 Recommended printer settings (strip printer)

Printer

Model	Char Set	Auto baudrate	Balance feature
P-2x	IBM/DOS	Off	Disable
P-52RUE	IBM/DOS ¹⁾	_	_
P-56RUE / P-58RUE	IBM/DOS ¹⁾ ²⁾	_	_

65

Instrument printer

Model	Baudrate	Bit / Parity	Stop Bits	Handshake
P-2x	9600	8/NO	1	Xon/Xoff
P-56RUE / P-58RUE	9600	8/NO	1	Xon/Xoff
	2)	2)	2)	2)
P-52RUE	9600	8 / None	1	XOn/XOff
	2)	2)	2)	2)

¹⁾ Printer settings not available.

7.5.3.2.3 Export management

Navigation: Home > Settings > Application settings > Print and export management > Export management

If a USB flash drive is connected and activated, final results can be automatically exported. The export contains information to the instrument, time, date, method, result as well as the graphical curve.



Only the formats .csv and .pdf can be exported on the USB flash drive.

The following parameters can be defined:

Parameter	Description	Values
File name prefix	Defines the name of the new data file.	any
File type	Defines the type of the data file.	CSV* I PDF
Printout language	Defines the language for general printouts. Note No impact on test or adjustment printouts.	English Deutsch Italiano Français Español Português Polski Magyar Český Русский 日本語 中文 한국어/조선말
Export management	Export results automatically after measurement if activated.	Selection: □* I ☑
Export target	Defines the destination of data transfer.	Peripheral

Export target: Peripheral

Parameter	Description	Values
Location	Finds the location on the storage device.	Browse and select

^{*} Factory setting

7.5.4 Quality management

Navigation: Home > Settings > Quality management

In this menu you can define all settings related to the adjustment and Test of the instrument. Periodic adjustments or tests ensure accurate measurement results.

7.5.4.1 Test / adjustment settings

Navigation: Home > Settings > Quality management > Test / Adjustment settings

This section describes all the menu options available for defining the parameters relating to the adjustment and testing of your instrument.

For performing test and adjustments **see** [Test/Adjust ▶ Page 50].

²⁾ Value applicable for Brazilian Portugese.

7.5.4.1.1 **Equipment**

Navigation: Home > Settings > Quality management > Test / Adjustment settings > Equipment

This menu item defines the equipment which are required for the following tests:

- Temperature Adjustment
- Temperature test
- SmartCal test (optional)
- Weight adjustment external
- Weight test external



For performing temperature adjustment or temperature test, you need the optional temperature adjustment kit, see [Accessories and Spare Parts > Page 87].

Test weight 1 and Test weight 2

The following parameters can be defined:

Parameter	Description	Values
Test weight name	Defines a name that is in an easily recognizable form as an alternative to the ID. A name should be unique and unambiguous.	any
Nominal weight	Defines the nominal weight value of the test weight. For adjustments, the weight must be at least 30 g.	0.1000 g 201 g
Identification	Defines the identification (ID) of the weight. The ID of the weight is provided on the Weight Certificate. The ID may contain your company specific identification number. This ID is stored in the adjustment or test history and shows up on the printouts.	any

^{*} Factory setting

Temperature kit

The following parameters can be defined:

Parameter	Description	Values
Temperature kit name	Defines a name that is in an easily recognizable form as an alternative to the ID. A name should be unique and unambiguous.	any
Identification	Defines the identification (ID) of the Temperature kit. The ID is provided on the Certificate. The ID may contain your company specific identification number. This ID is stored in the adjustment or test history and shows up on the printouts.	any

External humidity sensor

The following parameters can be defined:

Parameter	Description	Values
	Defines the identification (ID) of the RHT sensor. This ID is stored in the test history and shows up on the printouts.	any

For performing SmartCal tests, you can use a testo RHT sensor for the automatical setting of temperature and humidity values.

7.5.4.1.2 Test settings

Navigation: Home > Settings > Quality management > Test / Adjustment settings > Test settings

Moisture Analyzer Software Description

Weight test external

The following parameters can be defined:

Parameter	Description	Values
Tolerance	Defines the measuring tolerance of the external weight.	0.0001 % - 0.1000 % (0.0010 %)*
Test weight	To choose a test weight.	Test weight 1 I Test weight 2

^{*} Factory setting

Temperature test

The following parameters can be defined:

Parameter	Description	Values
Temperature kit	Shows the name of the Temperature kit which was defined in the menu item Equipment .	none
Temperature 1	Enables or disables the lower test temperature.	OFF I ON (100 °C)
Tolerance	Defines the tolerance of the test temperature.	15 °C (3 °C)*
Temperature 2	Enables or disables the upper test temperature.	OFF I ON (160 °C)
Tolerance	Defines the tolerance of the test temperature.	15 °C (3 °C)*

^{*} Factory setting

SmartCal test

■ Note

There are 4 fixed SmartCal methods and 2 customizable SmartCal methods. All 6 methods are activated as default values but the methods can individually be enabled or disabled.

The following parameters can be defined:

Parameter	Description	Values
Lot number input	Defines the type of input for the Lot number. The Lot number can be found on the product label. Free input: Input is not required Guided input: Input is required.	ON* I OFF I Free input I Guided input *
Clear lot number after measurement	Enabled or disabled the function for clearing the Lot number after test measurement.	☑ (enabled) I ☐ (disabled)*
Use cSmartCal	Sets the control limits to the correct values for using of cSmartCal.	✓ (enabled) I ☐ (disabled)*
Temperature	Defines the test temperature. Note If temperature in customer SmartCal methods is set, the control limits will be reset to default (off).	70 °C 100 °C 130 °C 160 °C Custom temperature 40 °C230 °C (105 °C*)

Control limits	Defines the control limits. The values of the control limits are in dependence of the temperature setting. The recommended values are automatically set. The values can also be changed. The recommended values are (lower - upper limits): 70 °C: 3.2-4.4 %MCN (cSmartCal: 3.3 - 4.3 %MCN) 100 °C: 5.2-6.4 %MCN (cSmartCal: 5.3 - 6.3 %MCN) 130 °C: 7.4-8.8 %MCN (cSmartCal: 7.5 - 8.7 %MCN) 160 °C: 9.9-11.7 %MCN (cSmartCal: 10.0 - 11.6 %MCN) Note If temperature in customer SmartCal methods is set, the control limits will be reset to default (off).	values see explanation I OFF
----------------	--	---------------------------------

^{*} Factory setting

7.5.4.1.3 Adjustment settings

Navigation: Home > Settings > Quality management > Test / Adjustment settings > Adjustment settings Weight adjustment external

The following parameters can be defined:

Parameter	Description	Values
Test weight	Selects the predefined Test weight .	Test weight 1 I
	Test weight 1 / Test weight 2 = defined in the menu item	Test weight 2
	Equipment	

Temperature adjustment

The following parameters can be defined:

Parameter	Description	Values
Temperature kit	Shows the name of the Temperature kit which was defined in the	none
	menu item Equipment .	

7.5.4.1.4 Print settings

Navigation: Home > Settings > Quality management > Test / Adjustment settings > Print and output settings

This menu item allows you to set specific printer functions for test or adjustment printouts.

- The setting for printout language in this menu item has impact on test or adjustment printouts only. For setting the general printout language, see [Print management ▶ Page 64].
- Not all languages are compatible with any printer, and also depending on the settings of the printer, see Recommended printer settings (Strip printer).
- For installing and defining a printer, see [Print management ▶ Page 64].

The following parameters can be defined:

Parameter	Description	Values
Printout language	Defines the printout language for test or adjustment printouts.	English I Deutsch I Italiano I Français I Español I Português I Polski I Magyar I Český
Print results automatically after test/ adjustment	Activated or deactivated function.	☑ (enabled) I ☐ (disabled)*

69

^{*} Factory setting

7.5.5 System and data management

Navigation: Home > Settings > System and data management

This menu item allows you to export or import settings and methods to or from an external USB storage device, as well as to restore the system.

Note for using USB memory stick

- USB memory sticks up to 32 GB are supported.
- Recommended formatting: FAT32 (NTFS not supported).
- Max. cluster size: 32 KB.
- Max. file size: 32 MB.
- Make sure that the stick is fully inserted.
- Make sure that the write protection is deactivated if present.

7.5.5.1 Export / Import

Navigation: Home > Settings > System and data management > Export / Import

This menu item allows you to export or import settings and methods to or from an external storage device. These data can be transferred to other instruments of the same type and can be useful to configure multiple instruments with the same specifications.



- Software Version must be equal.
- Data transfer between different Moisture Analyzer models (e.g. from HX204 to HS153 or HC103) is limited. Features that are not supported by a model and data resulting from these features cannot be imported.
- Export / import of methods can also be performed from the menu Method Definition.

Export settings and methods to external storage devices...

The following parameters can be defined:

Parameter	Description	Values
File name	Defines the name of the new data file.	any
Location	Finds the location on the storage device.	Browse and select
Export selection	Defines the data for export.	AII* I ✓ User
	 ■ Note	management and user preferences I Instrument, application and QM settings I
	User management and user preferences contains: User management settings, user preferences	
	Instrument, application and QM settings contains: Instrument settings, application settings, quality management settings	✓ Methods (x1)

^{*} Factory setting

- An external storage device is connected e.g. memory stick.
- Settings are made.
- To start, tap [Export].To abort, tap [Cancel].

Import settings and methods from external storage device...

The following parameters can be defined:

Parameter	Description	Values
Import from file	Defines the data to be imported.	Browse and select
Location	Shows the location of the import file.	_

Import selection	Defines the data for import.	All* I ✓ User
		management and user
		preferences I
		✓ Instrument, appli-
		cation and QM settings I
		✓ Methods (x1)

^{*} Factory setting

- Import settings and methods from external storage devices is selected.
- An external storage device is connected e.g. memory stick.
- 1 Tap Browse and select
 - → Import from file appears.
- 2 Select Device and import file.
 - → Full menu of Import settings and methods from external storage device... appears.
- 3 Tap **Import selection** and select the data for import if needed.
- 4 To start, tap the **Import** button and follow the instructions in the message window.
 - To abort, tap [Cancel].

7.5.5.2 Back up / Restore

Navigation: Home > Settings > System and data management > Back up / Restore

This menu item allows you to create a recovery point of the system and to store them on an external storage device. We advise you to create a recovery point of the system regularly. So you are able to restore the system with all data in fault.



Note

- After restoring, the system is in backup status. That means, that newer data such as settings, methods or results are lost.
- Restore is only possible on the same instrument as the back up was carried out.

Backup

The following parameters can be defined:

Parameter	Description	Values
File name	Defines the name of the new data file.	any
Location	Finds the location on the storage device.	Browse and select

- An external storage device is connected e.g. memory stick.
- Back up your system on external storage device... is selected.
- 1 Define the name of the new data file.
- 2 Define the location on the storage device.
- 3 To confirm, tap **OK**.
- 4 To start, tap **Backup**. To abort, tap **Cancel**.

Restore

The following parameters can be defined:

Parameter	Description	Values
Restore from file	Finds the location on the storage device and selects the file for	Browse and select
	restoring.	

71

- Restore your system from selected back up... is selected.
- An external storage device is connected e.g. memory stick.
- 1 Tap Restore from file... (Browse and select).

Moisture Analyzer Software Description

- → Storage device appears.
- 2 Select the file for restoring.
- 3 To start, tap **Restore**.
 - To abort, tap Cancel.
 - Confirmation window appears.
- 4 Confirm with **Restore**.
- The system will restore the data and then perform a restart. Wait until the system has completely started up again.

7.5.5.3 Export event logs

Navigation: Home > Settings > System and data management > Export event logs

This menu item allows you to export a log file (Zip format) with all relevant data to an external storage device. These data can be used for example, to support METTLER TOLEDO for troubleshooting.

The following parameters can be defined:

Parameter	Description	Values
File name	Defines the name of the new data file.	any
Export target	Defines the destination of data transfer.	Peripheral

Export target: Peripheral

Parameter	Description	Values
Location	Finds the location on the storage device.	Browse and select

^{*} Factory setting

7.5.5.4 Reset

Navigation: Home > Settings > System and data management > Reset

This menu item allows you to reset the Instrument to factory settings. Date, time, weight and temperature adjustments/history are not affected.



NOTICE

Data loss after system reset.

After resetting, the instrument is in factory default state. All data such as settings, methods or results is lost.

Save the data on an external storage device before you reset the instrument, see
 [Export / Import ▶ Page 70] and [Back up / Restore ▶ Page 71]

The following parameters can be defined:

Parameter	Description	Values
Reset	Resets the instrument to factory settings.	none

7.5.5.5 Update

72

Navigation: Home > Settings > System and data management > Update

The Update menu item is only used during the software update process, **see** [Updating the software ▶ Page 77].

Note

This function is not available on approved instruments.

Software Description Moisture Analyzer

7.5.6 Help and tutorials

Navigation: Home > Settings > Help and tutorials > Instrument tutorial > Next

The function **Instrument tutorial** automatically appeared when the function **Setup wizard** had been finished. The tutorial helps you, to define the basic functionalities of your Moisture Analyzer.

The application **Instrument tutorial** can be started directly at the end of the application **Setup wizard** or from the **Home** screen via the section **Settings**. The sections shown can be defined using the tutorial.

Moisture Analyzer Software Description

73

8 Maintenance

To guarantee the functionality of the Moisture Analyzer and the accuracy of the results, a number of maintenance actions must be performed by the user.

8.1 Maintenance tasks

Maintenance action	Recommended interval	Remarks
Cleaning	Depending on the degree of pollution or your internal regulations clean the instrument:	see chapter "Cleaning"
	After every use	
	After change of sample	
Performing routine tests	After cleaning	see chapter "Test"
(Weight test, temperature test, SmartCal test)	After a software update	
Performing adjustments	After changing the location	see chapter "Adjustments"
(weight adjustment, temperature adjustment)	If a test indicates that an adjustment is needed	
Replacing the dust filter (if used)	Depending on the degree of pollution	see chapter "Replacing the dust filter"
Replacing the power line fuse	If the fuse is blown	see chapter "Replacing the power line fuse"
Updating the software	If recommended by METTLER TOLEDO	see chapter "Updating the software"

8.2 Cleaning



MARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to injury and death. If the instrument cannot be shut down in an emergency situation, people can be injured or the instrument can be damaged.

 Disconnect the instrument from the power supply before cleaning or performing other maintenance work.



CAUTION

Burns due to hot surfaces

The interior parts of the heating module as well as the parts in the sample chamber can reach temperatures that can cause injuries if touched.

 Wait until the heating module has cooled down completely before performing any maintenance task.

Maintenance Moisture Analyzer

NOTICE

Damage to the instrument due to inappropriate cleaning agents

Inappropriate cleaning agents can damage the housing. If liquids enter the housing they can damage the instrument.

- 1 Make sure the cleaning agent is compatible with the material of the part you want to clean.
- 2 Make sure that no liquid enters the interior of the instrument. Do not spray any liquids and wipe off any spills immediately.
- 3 Use a lint-free cloth for cleaning.
- 4 Never open the housing of the instrument.

Protective equipment:


- Gloves
- Goggles

See also

Temperature adjustment ▶ Page 56

8.2.1 Sample chamber

- Sample chamber is open.
- 1 Remove the sample pan holder (1), the sample pan handler (2) and the draft shield element (3) for cleaning.
- 2 Carefully remove any deposits from the black temperature sensor (4).

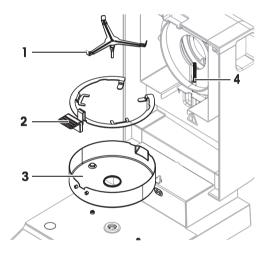

75

8.2.2 Heating module



↑ CAUTION

Burns due to hot surfaces


The round halogen lamp can reach temperatures that can cause injuries if touched.

- 1 Do not remove the halogen lamp.
- 2 Wait until the heating module has cooled down completely before cleaning.
- 3 Remove any splashes, deposits, or spots from the halogen lamp with a mild organic solvent such as ethanol.

Moisture Analyzer Maintenance

Overview

- 1 Reflector ring
- 2 Protective glass
- 3 Halogen lamp

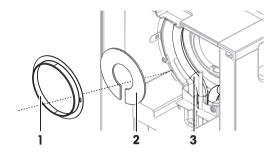


Removing reflector ring

Note

When the reflector ring is removed, the loose protective glass might fall out. Be careful when pulling out the reflector ring.

- 1 Sample chamber is open.
- 2 Unlock the reflector ring by pushing the anchoring clips on the side of the reflector ring inwards.
- 3 Carefully pull out the reflector ring.

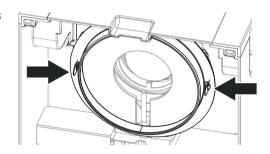


Removing protective glass

- Reflector ring is removed.
- Carefully pull out the protective glass.

Reassembling after cleaning

- All parts are cleaned.
- 1 Insert the protective glass.
- 2 Set the protective glass into the housing and adjust the position until it sits correctly.
- 3 Insert the reflector ring and push the anchoring clips into the designated openings.
- 4 Close the heating module.

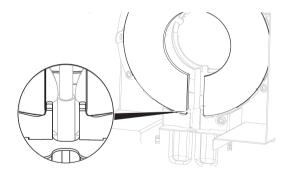

8.2.3 Fan grill

76

The air inlet of the fan is located at the rear of the instrument and its exterior should be cleaned from time to time to free it from any dust deposits.

8.2.4 Putting into operation after cleaning

- 1 Reassemble the Moisture Analyzer.
- 2 Press **(**) to switch on the Moisture Analyzer.
- 3 Warm up the Moisture Analyzer. Wait 1h for the acclimatization, before starting the tests.
- 4 Check the level status, level the Moisture Analyzer if necessary.


Maintenance Moisture Analyzer

- 5 METTLER TOLEDO recommends to perform weight and temperature tests and, if necessary, adjustments after cleaning the Moisture Analyzer.
- → The Moisture Analyzer has been put into operation and is ready to use.

8.3 Replacing the dust filter

If you use a dust filter for your Moisture Analyzer, check the filter pollution at regular intervals. For replacement filters **see** Accessories and Spare Parts.

- 1 Open the filter lid.
- 2 Replace the dust filter.

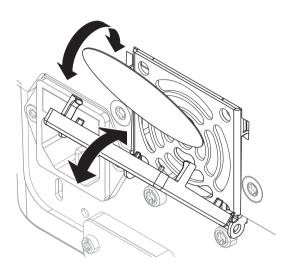

77

8.4 Replacing the power line fuse



MARNING

Danger of death or serious injury due to electric shock!

Risk of electrical shock or damage of the instrument when changing fuses.

- 1 Do not use a fuse of a different type or rated value.
- 2 The short-circuiting (bridging) of fuses is not allowed.

The power line fuse is located on the back of the drying unit. To change the fuse, proceed as follows:

- 1 Pull out the power cord.
- 2 Slide out the fuse holder (1) with a suitable tool such as a screwdriver.
- 3 Remove the fuse (3) and check its condition.
- 4 If the fuse is blown, replace the fuse with one of the same type and the same rated value.

 Note

A spare fuse is located in the holder (2).

8.5 Updating the software

METTLER TOLEDO is constantly improving its instrument software to make the handling of its instruments more comfortable and smoother. The current software versions can be downloaded online and installed by the user to benefit from the improvements quickly and easily.

The firmware made available on the Internet has been developed and tested by Mettler-Toledo GmbH using processes that meet the guidelines of ISO 9001. Mettler-Toledo GmbH does not, however, accept liability for consequences that might arise from using the firmware.

Moisture Analyzer Maintenance

NOTICE

Data loss during software update.

During the update of the software, data such as settings, methods or results can be lost. This is also the case if the update fails.

 Before starting a software update, create a backup file of the current system on an external storage device.

Search for software

► http://www.mt.com/moisture-software

Update procedure

- Backup of the current data is created, see [Back up / Restore ▶ Page 71]
- Recent software update is downloaded and stored on an external device, e.g., USB stick or SD card.
- Administrator rights are active.
- 1 Insert device with software for update into its respective port.
- 2 Wait until the external device is recognized by the instrument.
- 3 Select Home > Settings > System and data management > Update. If Update is not visible check if administrator rights are active. If you have an approved version, contact your support representative.
 - → Update screen appears.
- 4 Select the device on which the software is stored.
- 5 Confirm with [**Update**] and wait until the software update is completely finished. Do not remove the device before the instrument has completely restarted.
 - → The instrument reboots.
 - → A message appears: Operating system is up to date. Update instrument software?...
- 6 Confirm with [Update].
- 7 Wait until the system returns to the home screen.

Software downgrade

It is not recommended to downgrade the software. However, in case of a setup that requires a software downgrade in order to work, please observe the following safety instructions.



NOTICE

Data loss during software downgrade

During a downgrade of the software all device data such as drying methods, results, SmartCal history, and settings will be removed.

Only let a service technician perform a software downgrade.

Maintenance Moisture Analyzer

9 Troubleshooting

Possible errors with their cause and remedy are described in the following chapter. If there are errors that cannot be corrected through these instructions, contact METTLER TOLEDO.

9.1 Error messages

Error message	Possible cause	Diagnostic	Remedy
Underload	Underload - the sample pan holder is missing.	Check.	Insert the sample pan holder. If needed, reboot the system by discon- necting and reconnecting the power supply.
Overload	Overload - the weight on the sample pan exceeds the weighing capacity of the balance.	_	Reduce the weight of the sample.
Blinking zeros	The balance is out of the zero range. When the instrument was switched on or upon zeroing, one or more limits were exceeded. The usual reason for this is when there is a weight on the weighing pan when the balance is switched on.	Check if a weight is on the sample pan.	Remove the weight and retry.
"Unhandled exception caught in main thread."	Startup problem on terminal.	_	Download and install the appropriate software. See [Updating the software ▶ Page 77].
"Sensor connection lost!"	More than 5000 measurements have been stored on the memory.	With software V2.11 and earlier, this caused the sensor to lose connection.	Update the software to V2.12 or later, see [Updating the software ▶ Page 77]

9.2 Error symptoms

Error symptom	Possible cause	Diagnostic	Remedy
Moisture analyzer cannot be switched on.	The power supply cable is not connected.No power supply voltage in the mains.	Check.	Connect the power supply cables or reestablish the mains supply voltage.
	Faulty fuse.	Check.	Replace the fuse. The fuse is located on the back of the instrument besides the power plug (find a spare fuse there).
Keys and buttons on the terminal do not function.	Software bug.	_	Restart the software by disconnecting and reconnecting the power plug.
Weight out of range. During adjustment.	Either no weight piece or the wrong weight piece has been placed on the	Check if the correct weight piece is placed.	Use the correct weight piece and retry.

79

Moisture Analyzer Troubleshooting

Error symptom	Possible cause	Diagnostic	Remedy
	sample pan for adjustment. This message is also displayed if you do not remove the weight piece when you are prompted.		
Drift The measuring values on the display slowly drift up/down.	The air current which cools the Load cell is blocked (filter, grating). The rising temperature influences the	Check (find the air filter and the grating at the backside of the moisture analyzer).	Replace filter (if available) and clean ventilation grating.
Unstable weight indication.	measurement. The sample pan holder is not placed properly.	Check.	Make sure the sample pan holder is correctly placed.
The measuring values do not stabilize but go up/			Some parts are touching the sample pan.
down.			Volatile sample substance (weight of the sample changes quickly).
The measurement takes too long.	An unsuitable switch-off criterion was set.	_	Choose a suitable switch-off criterion.
	Sample substance tends to form a skin, when heated.		If you use samples which tend to form a skin that hinders evaporation, perform the measurement at a higher temperature.
	An excessive amount of sample cause slow drying.	_	Enlarge the surface of the sample substance, e.g by crushing or grinding.
	Liquids take longer to dry.	_	For sample liquids, use absorbent glass fibre filters.
			Use absorbent glass filter for liquids.
			Enlarge the surface of the sample, e.g. by crushing or grinding.
Measurement results are not repeatable.	Unstable environment/ location of the instrument.	_	Choose a suitable location, see [Selecting the location ▶ Page 12].
			The drying time is too short for the "Timed switch-off" criterion.
	The sample substance boils and the splashed drops continuously change the weight.	_	Lower the drying temperature.
			The samples are not comparable.
	The granulation of the sample is not homogenous or too large.	_	Use samples with a homogenous granulation.

Troubleshooting Moisture Analyzer

Error symptom	Possible cause	Diagnostic	Remedy
	Insufficient heating power because the protective glass of the halogen radiator is dirty.	Check if the protective glass is dirty.	• Clean the protective glass, see [Cleaning • Page 74].
	The temperature sensor is contaminated/dirty.	Check if the temperature sensor is dirty.	Clean the temperature sensor, see [Cleaning ▶ Page 74].
	The sample substance does not become completely dry due to uneven distribution in the sample pan.		Evenly spread the sample substance in the pan and retry.
The difference between target and actual temperature is out of tolerance. When using the temperature adjustment kit.	The temperature adjustment kit is not adjusted or defective.		 Try another temperature adjustment kit. Adjust the temperature adjustment kit (contact your METTLER TOLEDO service representative)
The adjustment temperature is lower than expected (e.g. 90°C instead of 100°C).	The draft shield is not installed.	Check.	Install the draft shield.
The SmartCal test failed.	The test prerequisites are not met correctly.	_	Make sure to meet the test prerequisites, see [SmartCal test > Page 53].
	The entered room temperature and/or the	Check if the values are set correctly.	Enter the correct values and retry.
	relative humidity are set wrong.		The test substance does not become completely dry due to uneven distribution in the sample pan.
	The test substance does not become completely dry due to uneven distribution in the sample pan.	_	Evenly spread the test substance in the pan and retry.
	The instrument is out of adjustments.	_	Perform:
			Temperature test, see Temperature test.
			2. Weight tests, see Tests
			3. Repeat the SmartCal test with a new test substance.

Moisture Analyzer Troubleshooting

81

10 Technical Data

10.1 General data

Power supply

110 V AC Version 100 V–120 V, 50/60 Hz, 4 A 230 V AC Version 200 V–240 V, 50/60 Hz, 2 A

Voltage fluctuations -15%+10%

Power load max. 450 W during drying process

Power line fuse $115 \text{ V: } 5 \times 20 \text{ mm, } \text{F6.3 AL } 250 \text{ V } \text{(6.3 A, fast-acting, low-$

breaking capacity)

230 V: 5×20 mm, F2.5 AL 250 V (2.5 A, fast-acting, low-

breaking capacity)

Protection and standards

Overvoltage category II
Degree of pollution 2

Standards for safety and EMC see Declaration of Conformity (part of standard equipment)

Range of application for use in dry interior rooms

Environmental conditions

Height above sea level up to 4000 m

Ambient temperature range Operation: +10 °C to 30 °C

(operability guaranteed 5 °C to 40 °C)

Relative air humidity max. 80% up to 31 °C, linearly decreasing to 50% at 40 °C

20% - 80% and non-condensing conditions

Warm-up time At least 60 minutes after connecting the instrument to the power

supply; when switched on from standby, the instrument is ready

for operation immediately.

Materials

Drying Unit

Housing PBT, PBX45A (UL 94 V-0) Inspection window grill PPS, A504X90 (U L94 V-0)

Protective glass
Halogen lamp
Quartz glass

Reflector bracket PPS A504X90 (UL 94 V-0)

Draft shield, interior bottom plate Stainless steel, X2CrNiMo17-12-2 (1.4404)

Technical Data Moisture Analyzer

10.2 Model-specific data

Drying unit

Heating module Halogen ring-shaped radiator

Temperature range 40–230 °C
Temperature step 1 °C

Temperature programs standard, rapid Drying time 480 minutes

Balance

Maximum capacity 101 g
Readability 1 mg
Minimum sample weight 0.5 g
Weighing technology Strain gauge

Adjustment External weight (100 g, accessory)

Moisture Content

Readability 0.01%
Repeatability (sd) with 2 g sample 0.10%
Repeatability (sd) with 10 g sample 0.015%

Interfaces

1 × RS232C (for production only)1 × RS232C (9-pin socket)

2 × USB Host (Type A socket), USB 1.1

Supported memory sticks up to 32 GB

1 × USB Device (Type B socket), USB 1.1

83

Data

Export data format of results
• CSV (Comma-separated values)

PDF

Hardware

Drying unit

Sample chamber opening/closing manual

Leveling 2 leveling screws

Sample pan ø 90 mm Maximum height of sample 25 mm

Thermal overload protection Bimetallic-element switch in heating module

Dimensions (w \times h \times d) 192 \times 169/370 \times 385

Weight, ready to measure 5.1 kg


Display WVGA color display, resistive touch-sensitive

Moisture Analyzer Technical Data

10.3 Dimensions

84

(all dimensions in mm)

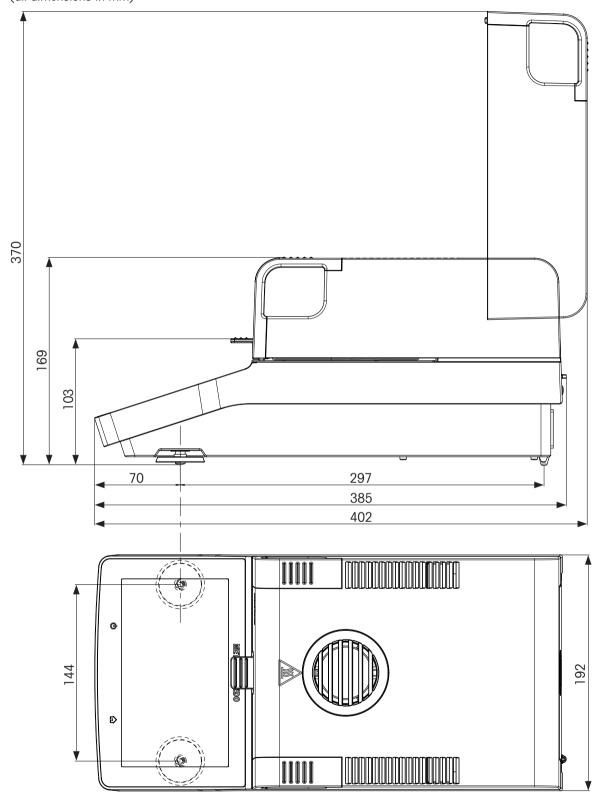

Technical Data Moisture Analyzer

10.4 Interfaces specification

10.4.1 RS232C

Schematic	Item	Specification
	Interface type	Voltage interface according to EIA RS-232C/ DIN66020 CCITT V24/V.28)
DATA	Max. cable length	15 m
RXD IN TXD OUT	Signal level	Outputs: +5 V +15 V (RL = 3-7 k Ω) -5 V15 V (RL = 3-7 k Ω) Inputs: +3 V +25 V -3 V25 V
5	Connector	Sub-D, 9-pole, female
	Operating mode	Full duplex
90 0 0 6 HAND	Transmission mode	Bit-serial, asynchronous
CTS SHA KE	Transmission code	ASCII
RTS	Baud rates	see setting options
OUT	Bits/parity	see setting options
	Stop bits	see setting options
	Handshake	None, XON/XOFF, RTS/CTS (software selectable)

10.4.2 USB host

Schematic	Item	Specificati	on	
	Standard	In conform 1.0/1.1	ity with USB Specification Revision	
	Speed	Full speed	Full speed 12 Mbps (requires shielded cable)	
] [Power usage	Max. 500	mA	
	Connector	Type A		
1 2 3 4	Pin assignment	1	VBUS (+5 V DC)	
		2	D- (Data -)	
		3	D+ (Data +)	
		4	GND (Ground)	
		Shell	Shield	

Moisture Analyzer Technical Data

10.4.3 USB device

Schematic	Item	Specification
2 1	Standard	In conformity with USB specifications revision 2.0
	Speed	Full-speed 12 Mbps (requires shielded cable)
	Function	CDC (Communication Device Class) serial port emulation
3 4	Power usage	Suspended device: Max 10 mA
	Connector	Type B
1 VBUS (+5 VDC)		
2 D- (Data -)		
3 D+ (Data +)		
4 GND (Ground)		
Shield Shield		

Technical Data Moisture Analyzer


11 Accessories and Spare Parts

11.1 Accessories

	Description	Order no.
Sample Handling		
	Aluminum sample pans, set of 80 pcs	00013865
	Glass fiber filter (for Liquids), set of 100 pcs	00214464
	Extra strong aluminum sample pans, set of 80 pcs	11113863
	Reusable steel sample pans, height 6 mm, set of 3 pcs	00214462
	Textile cage for bulky samples (HA-cage), 1 pcs	00214695
	Down holder, prevents slipping of the sample during measurement, 1 pcs	00214758
Quality Management		
	Certified adjustment weight, 100 g (F1)	30406400



OIML / ASTM weights (with calibration certificate) see www.mt.com/weights

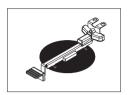

Certified temperature kit 30020851



SmartCal Moisture Analyzer Test Substance

cSmartCal, certified, 12 tests	30005793
cSmartCal, certified, 24 tests	30005791
SmartCal, 12 tests	30005792
SmartCal, 24 tests	30005790
StarterPac cSmartCal	30005918
StarterPac SmartCal	30005917

Printers



RS-P25/01 (EMEA) printer with RS232C connection to 11124300 instrument

Paper roll (length: 20 m), set of 5 pcs 00072456

Paper roll (length: 13 m), self-adhesive, set of 11600388
3 pcs

00065975

11600388

30237290

30094673

Ribbon cartridge, black, set of 2 pcs



USB-P25/01 (EMEA) printer with USB connection to instrument
Paper roll (length: 20 m), set of 5 pcs
11124301
00072456

Paper roll (length: 13 m), self-adhesive, set of 3 pcs

Ribbon cartridge, black, set of 2 pcs 00065975



P-52RUE dot matrix printer RS232C, USB and Ethernet connections, simple print-outs

Paper roll (length: 20 m), set of 5 pcs 00072456

Paper roll (length: 13 m), self-adhesive, set of 11600388

3 pcs

Ribbon cartridge, black, set of 2 pcs 00065975



P-56RUE thermal printer with RS232C, USB and Ethernet connections, simple print-outs, date and time

Paper roll, white (length: 27 m), set of 10 pcs 30094723 Paper roll, white, self-adhesive (length: 13 m), set 30094724

of 10 pcs

Accessories and Spare Parts

Moisture Analyzer



P-58RUE thermal printer with RS232C, USB and Ethernet connections, simple print-outs, date and time, label printing, balance applications, e.g., statistics, formulation, totaling

> Paper roll, white (length: 27 m), set of 10 pcs 30094723 Paper roll, white, self-adhesive (length: 13 m), set 30094724 of 10 pcs

> Paper roll, white, self-adhesive labels (550 labels), 30094725 set of 6 pcs

Dimension of the label 56×18 mm

Anti-theft devices



Anti-theft cable with lock

11600361

30094674

Cables for RS232C interfaces



RS9 – RS9 (m/f): connection cable for PC, length = 1 m

11101051

Barcode readers



Corded USB barcode reader

30417466


Cables for USB interface



USB (A -B) connection cable, length = 1 m


30241476

Miscellaneous

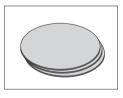


Dust filter, set of 50 pcs

11113883

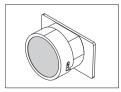


HC dust filter housing

30216118

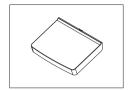

Software

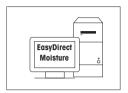


EasyDirect Moisture (simple data collection)

30522676

11.2 Spare parts

90

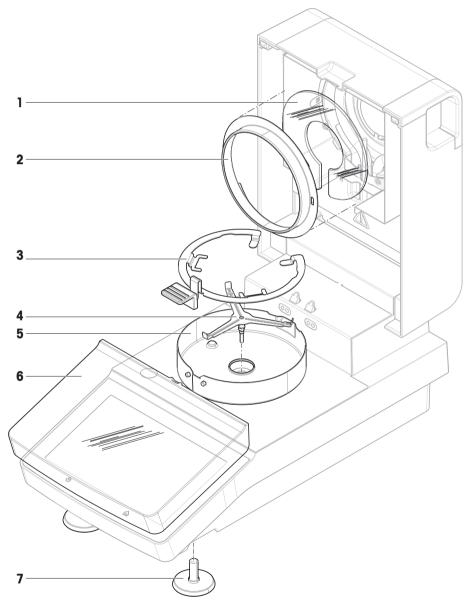

	Order no.	Designation	Remarks
1	30216116	Protective glass	_

Accessories and Spare Parts

Moisture Analyzer

	Order no.	Designation	Remarks
2	30216117	Reflector ring	_
3	30216113	Sample pan handler	_
4	214642	Sample pan holder	_
5	30216114	Draft shield	_
6	30216115	Protective cover display	_
7	30104835	Pair of leveling feet	_

12 Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.



Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties, the content of this regulation must also be related.

Disposal Moisture Analyzer

Index

Accessories 37, 87 EasyDirect Moisture 48 Adding comments 43 Edit Adjustment 16, 22 Method 40 Balonce 55 Entering Heding module 55 Numerical values 22 Settlings 66 Text and numbers 23 Settlings 66 Export 72 Anti-theff 15 Export 72 Asterisk 34, 36 Export 72 Asterisk 34, 36 Export 72 Automatic sample chamber 37 Resull during and methods 73 B Forth Forth 72 Asterisk 34, 36 Export 48, 8 Brighthess 60 F 6 C C Forth 6 6 C C Forth 6 76 76 77 76 77 76 77 76 77 76 77 72 <	A		E	
Adjushment 16, 22 Method 40 Balance 55 Entering 24 Heatling module 55 Numerical values 24 Settings 66 Text and numbers 22 Settings 66 Text and numbers 22 Weight external 55 Event logs 77 Anti-theft 15 Export 72 Asterisk 34, 36 Export 48, 83 B Settings and methods 77 Back up 71 Result 48, 83 Settings and methods 77 External weight adjustment 68 C C Factor 36 Cortified RHT sensor 54 Filter 44 Certified RHT sensor 54 Filter 47 Chord 45 Fire factor 36 Cortified RHT sensor 43 Filter 47 Chord 48 Filter 42 Cortified RHT sensor 48 <t< th=""><th>Accessories</th><th>37, 87</th><th>EasyDirect Moisture</th><th>49</th></t<>	Accessories	37, 87	EasyDirect Moisture	49
Bolance 55 Entering 24	Adding comments	43	Edit	
Heating module 55	Adjustment	16, 22	Method	40
Seffings 66 Text and numbers 23 Touch screen 63 environmental condition 12 Weight external 55 Event logs 72 Anti-theff 15 Export 72 Asterisk 34, 36 Export 48, 83 Automatic sample chamber 37 Result 48, 83 Back up 71 External weight adjustment 68 Brightness 60 F Foctor 36 Brightness 60 F Foctor 36 Certified RHT sensor 54 Filter 48 Cortified RHT sensor 54 Filter 48 Comments to measurement 43 Free factor 36 Comments to measurement 43 General method properties 37 Coty 83 General method properties 37 CSV 83 General method properties 37 CSV format 48 Graphics view 48 Date 25	Balance	55	Entering	
Touch screen	Heating module	55	Numerical values	24
Weight external 55 Export 72 Anti-theff 15 Export 72 Asterisk 34, 36 Export 72 Asterisk 34, 36 Export 48, 83 Automatic sample chamber 37 Result 48, 83 Back up 71 External weight adjustment 68 Back up 71 Factor 36 Back up 71 Factor 36 Colculated values 34 Factor 36 Certified RHT sensor 54 Filter 47 Chotral 45 Free factor 36 Comments to measurement 43 Free factor 36 Comments to measurement 43 Free factor 36 Covy 83 General settings 22 CSV 83 General settings 22 CSV format 83 H Data management 70 History 56 Data management 70 <t< td=""><td>Settings</td><td>66</td><td>Text and numbers</td><td>23</td></t<>	Settings	66	Text and numbers	23
Anti-theff 15 Export 72 Asterisk 34, 36 Export Automatic sample chamber 37 Result 48, 83 B Settings and methods 77 Back up 71 External weight adjustment 68 Brightness 60 F C Factor 36 Colladidated values 34 Filter Certified RHT sensor 54 Filter 45 Chart 45 Free factor 36 Comments to measurement 43 Free factor 36 Compt G G Method 41 General method properties 37 CSV 83 General settings 22 SV format 48 Graphics panel 22 CSV format 48 History 56 Data 83 Hence screen 18, 20, 22 Define Host settings 61 Method 39, 40 Hence screen	Touch screen	63	environmental condition	12
Asterisk 34, 36 Export Automoffic sample chamber 37 Result sellings and methods 37 BC Seltings and methods 37 Brightness 60 F C Factor 36 Calcidulated values 34 Filter 45 Certified RHT sensor 54 Filter 45 Chord Free factor 36 Comments to measurement 43 Free factor 36 Comments to measurement 43 Free factor 36 Copy 6 6 Welthod 41 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics view 48 Data 83 H H Data 83 H H Date 25 Home screen 18, 20, 22 Define 40 Home screen 18, 20, 22 Method 39, 40 Identi	Weight external	55	Event logs	72
Automatic sample chamber 37 Result 48, 85 B Settings and methods 70 Back up 71 External weight adjustment 68 Brightness 60 F C Factor 36 Calciulated values 34 Filter 45 Certified RHT sensor 54 Filter 45 Chard 45 Free factor 36 Comments to measurement 43 Free factor 36 Comy General method properties 37 36 Coty 83 General settings 32 CSV format 48 Graphics view 45 Data management 70 History 56 Data management 70 History 56 Define 45 Horrescreen 18, 20, 22 Define 45 Horrescreen 18, 20, 22 Define 45 Horrescreen 18, 20, 22 Define 45 Horrescreen <t< td=""><td>Anti-theft</td><td>15</td><td>Export</td><td>72</td></t<>	Anti-theft	15	Export	72
Book up 71 Brightness Settings and methods 70 External weight adjustment 70 External weight adjustment 70 External weight adjustment 66 External weight adjustment 70 External weight adjustment 66 External weight adjustment 70 External weight adjustment 66 External veight adjustment 67 External veight adjustment 68	Asterisk	34, 36	Export	
Back up 71 External weight adjustment 68 Brightness 60 F C Factor 36 Calculated values 34 Filter 45 Certified RHT sensor 54 Filter 45 Chart 45 Free factor 36 Comments to measurement 43 Free factor 37 Cony G G Method 41 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics ponel 25 CSV format 48 Graphics ponel 25 Data 83 H History 58 Date 25 Home screen 18, 20, 22 22 Define 40st settlings 61 61 Method 39, 40 Identifications 63 Display in mode 34 Identifications ponel 23 Display mode 34 Identificati	Automatic sample chamber	37	Result	48, 83
External weight adjustment Section F	В		Settings and methods	70
Brightness 60 F C Factor 36 Calculated values 34 Filter 45 Certified RHT sensor 54 Filter 45 Chart 45 Firee factor 36 Comments to measurement 43 Free factor 36 Comments to measurement 43 Fuse 77, 82 Convention 5 G Copy 6 G G Method 41 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics view 45 Data management 70 History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 Identifications 63 Delete see Remove 41 Identifications panel 23 Display longuage 59 Settings and methods 70		71	External weight adjustment	69
C Cactor Factor 38 Calculated values 34 Filter 45 Certified RHT sensor 54 Filter 45 Chard 45 Free factor 36 Comments to measurement 43 Free factor 36 Convention 5 Free factor 36 Copy 6 6 Method 41 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics panel 25 CSV format 83 H H Data 83 History 56 Data 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 Interface 63 Delete 1 Identifications panel 23 See Remove 41 Identifications panel 23 Disposal 92 Settings and methods 70	•		F	
Fan 76, 77 Calculated values 34 Filter 48 Certified RHT sensor 54 Filter 48 Certified RHT sensor 54 Filter 48 Comments to measurement 43 Filter 77 Comments to measurement 43 Filter 48 Free factor 38 Free factor 38 Fuse 77 Suse	Brigniness	60		20
Calculated values 34 Filter 45 Certified RHT sensor 54 Firmware update 77 Chard 45 Free factor 36 Comments to measurement 43 Free factor 36 Convention 5 6 Copy 6 6 Method 41 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics panel 23 Orata 83 H History 58 Date 25 Home screen 18, 20, 22 Define 45 Host settings 61 Method 39, 40 Intensions 63 Define 41 Identifications 63 Method 39, 40 Intensions 63 Display longuage 59 Identifications 63 Display mode 34 Identifications 63 Drying 50 Settings and met	C			
Certified RHT sensor 54 Chart Firmware update 77 Chart 45 Free factor 36 Comments to measurement 43 Fuse 77, 82 Convention 5 6 Copy 6 6 Method 41 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics panel 23 Data 83 H H Data management 70 History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 I Identifications 63 Method 39, 40 I Identifications 63 Display language 59 Settings and methods 70 Display language 59 Settings and methods 70 Display mode 34 Settings and methods 70 Display program Interface 83 Rapid 31 Introduction 83 Total method 84 Interface	Calculated values	34		•
Chart 45 Free factor 36 Comments to measurement 43 Free factor 36 convention 5 G Copy G 37 Method 41 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics panel 23 Dotal 83 H Data 83 H Data 83 H Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 I Delete See Remove 1 Dimensions 84 Identifications 63 Display Inguage 59 Settings and methods 70 Display mode 34 Install 12 Drying Interface 83 Drying program Interface 83 Introduction 84 Introduct	Certified RHT sensor	54		
Convention 5 Fuse 77, 82 Convention 5 G Copy General method properties 37 Method 41 General settings 22 CSV 83 General settings 22 CSV format 48 Graphics panel 25 Dotata 83 H History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 I Delete Identifications 63 see Remove 41 Identifications panel 23 Display language 59 Settings and methods 70 Display mode 34 Install 12 Drying install 12 Temperature 31, 32 Interface 83 Drying program Interface Interface 83 Interface Interface 10 Interface Interface 1	Chart	45	•	
Convention Copy Method A11 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics panel 23 Data Bata B	Comments to measurement	43		
Method 41 General method properties 37 CSV 83 General settings 22 CSV format 48 Graphics panel 23 Data 83 H Data management 70 History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 I Delete Identifications 63 see Remove 41 Identifications panel 23 Display language 59 Settings and methods 70 Display mode 34 Settings and methods 70 Disposal 92 install site 12 Drying site Interface 83 Drying program Interim results 64 Interim results Introduction 84 Introduction 85	convention	5		11, 02
CSV 83 General settings 22 CSV format 48 Graphics panel 23 D Graphics view 45 Data 83 H Data management 70 History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 I Delete Identifications 63 see Remove 41 Identifications panel 23 Display language 59 Settings and methods 70 Display mode 34 Install 31 Drying 11 12 Temperature 31, 32 Interface 83 Drying program Interface 83 Rapid 31 J	Сору		G	
CSV format 48 Graphics panel 23 Data 83 H Date management 70 History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 Identifications 63 Dielete Identifications panel 23 Dimensions 84 Import 23 Display language 59 Settings and methods install 70 Disposal 32 Interface 83 Drying Interface 83 Drying program Interface 83 Rapid 31 J	Method	41	General method properties	37
Data 83 H Date management 70 History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 I Dielete Identifications 63 see Remove 41 Identifications panel 23 Display language 59 Settings and methods 70 Display mode 34 Settings and methods 70 Disposal 92 install site 12 Drying Interface 83 Drying program Interface 83 Rapid 31 Introduction 58 Standard 31 J	CSV	83	General settings	22
Data management 70 History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 Delete see Remove 41 Identifications panel 19, 23 Display language 59 Import 20, 23 Display mode 34 Settings and methods install 19, 20, 21 Drying Temperature 31, 32 Interface 19, 21 Drying program Rapid 31 Standard 31 J	CSV format	48	Graphics panel	23
Data management 70 History 58 Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 Delete see Remove 41 Identifications panel Import 23 Display language 59 Settings and methods install 25 Disposal 92 install 27 Drying Temperature 31, 32 Interface 83 Rapid Standard 31 J	D		Graphics view	45
Date Define Host settings 61 Method 39, 40 Delete see Remove Dimensions Display language Display mode Disposal Drying Temperature Drying program Rapid Rapid Standard 18, 20, 22 Home screen Host settings Host settings Host settings Host settings Identifications Host settings Identifications Host settings Identifications For identifications Host settings Identifications For identifica	Data	83	Н	
Date 25 Home screen 18, 20, 22 Define Host settings 61 Method 39, 40 Delete see Remove 41 Identifications panel Import 23 Display language 59 Settings and methods install 23 Drying Temperature 31, 32 Interface Interim results Introduction 54 Drying program Rapid 31 Standard 31 J	Data management	70	History	58
Define Method 39, 40 Delete see Remove Dimensions Display language Display mode Disposal Drying Temperature Drying regram Rapid Standard Method 39, 40 I Identifications Identifications panel Import Settings and methods install site Interface Interface Interface Interim results Introduction Introduction Introduction Introduction Introduction Introduction Introduct		25	Home screen	18, 20, 22
Method Delete See Remove See Remove Dimensions Display language Display mode Disposal Drying Temperature Drying rogram Rapid Standard 39, 40 Identifications Identifications panel Import Settings and methods install Interface Interface Interim results Introduction Identifications Identifications Interface Import Import Settings and methods Interface Interim results Interim results Introduction Introduction Identifications Identifications Import Import Import Interim results Interim results Introduction Introduction Introduction Identifications Import Import Import Interim results Interim results Introduction Introduction Introduction Identifications Import Import Import Interim results Interim results Interim results Introduction Introduction Introduction Interim results Introduction Interim results Introduction Introduction Interim results Introduction Interim results Introduction Interim results Introduction Interim results Introduction Introduction Interim results Interim results Introduction Interim results Introduction Interim results Interim results Introduction Interim results Interim r	Define		Host settings	
see Remove 41 Identifications 63 Dimensions 84 Identifications panel 23 Display language 59 Import Settings and methods 70 Display mode 92 install site 12 Drying Temperature 31, 32 Interface Interim results 64 Drying program Rapid 31 Standard 31 J	Method	39, 40	1	
Dimensions Display language Display mode Disposal Drying Temperature Drying program Rapid Standard All Identifications panel Import Settings and methods install site Interface Interface Interoduction Introduction Introduction Standard Introduction J	Delete			
Display language Display mode Disposal Drying Temperature Drying program Rapid Standard Display language Settings and methods install site Interface Interface Interim results Introduction Introduction Settings and methods Install site Interface Interim results Introduction Standard J	see Remove	41		
Display language Display mode Disposal Drying Temperature Drying program Rapid Standard Settings and methods install site Interface Interim results Introduction Standard Settings and methods install site Interface Interim results Introduction Standard J	Dimensions	84		23
Disposal 92 install Drying	Display language	59	·	
Drying Temperature 31, 32 Drying program Rapid Standard 31 J site 12 Interface Interim results Introduction 5 J	Display mode	34	_	/(
Temperature 31, 32 Interface 83 Drying program Rapid 31 Standard 31 J	Disposal	92		1.0
Drying program Rapid Standard 31, 32 Interim results Introduction 5 5 5 5 5 64	Drying			
Rapid 31 Introduction 5 Standard 31 J	Temperature	31, 32		
Standard 31 J	Drying program			
	Rapid	31	introduction	5
Dust filter 77 Journal 45	Standard	31	J	
	Dust filter	77	Journal	45

Moisture Analyzer Index 93

K		Power line fuse	//
Kensington slot	15	Power supply	82
Keyboard		Print management	64
text and numbers	23	Printer	0.7
Keyboard language	59	Settings	61
Keypad		Printout	
numerical values	24	Intermediate results	64
Keys on the terminal	18, 20	Language	64
L		Q	
Languages	59	QuickAccess	22
Printout	64	R	
Leveling	14	Rapid drying	31
Lists	25	Remove	
location	12	Method	41
Log file	72	Replace	
M		Dust filter	77
		Power line fuse	77
Maintenance	77	Reset	72
Dust filter	77	Restore	71
Power line fuse	77	Result	
Measurement	22	Export	48, 83
First measurement	26	Result and value handling	36
Simple measurement	26	Results	22, 45
Memory stick	18, 70	RHT sensor	54, 67
Method			2.7, 2.
Copy	41	S	
Definition	30	safety information	7
Edit	22, 40	Sample preparation	44
Export and import	70	Scope of delivery	12
Name	38	Settings	22
New	22, 39, 40	Adjustment	66
Properties	37	Date	25
Remove	41	Export and import	70
Renaming	38	Host	61
MethodWizard	38	Peripheral	61
MT-SICS	61	Temperature Test	68
N		Time	25
New method	39, 40	Weight adjustment	69
Numbers	23	Weight test	68
Numerical values	24	Shortcuts	21, 22, 23, 44
P		Show results	22
	00	SmartCal	F0
Parameter panel	23	General	53
Password	60	Test cettings	54
PDF	83	Test settings	68
Peripheral settings	61	SOC	32

Index Moisture Analyzer

Software update	77
Sound	60
Standard drying	31
Standard equipment	12
Standby	18, 20
Start measurement	22
Start weight	35, 44
Starting mode	37
Status messages	21
Stop drying	43
Switch button	24
Switch ON/OFF	18, 20
Switch-off criterion	32
symbol	5
warning	7
System management	70
T	
Tables	25
Temperatur	
Adjustment settings	69
Temperature	
Switch-off criterion	32
Test	52
Test settings	68
Test	22
Balance	51
Heating device	51
Temperature	52
Weight external	53, 55
Tests	
General	51
Text	23
Time	25
Touch screen adjustment	63
Troubleshooting	72
U	
Unpacking	12
USB Driver	
Installation	17
USB Memory stick	
see memory stick	18, 70
User	
Home screen	22
Interface	20
Shortcuts	22

V	
Value panel	23
W	
warning symbol	7
Weighing-in aid	36, 44
Weight	
Adjustment external	55
Test external	53, 55
Weight adjustment	
Settings	69
Weight test	
external	53
Settings	68
Work screen	23
Workflow handling	37

Moisture Analyzer Index 95



Good Weighing Practice[™]

GWP® is the global weighing standard, ensuring consistent accuracy of weighing processes, applicable to all equipment from any manufacturer It helps to:

- Choose the appropriate balance or scale
- Calibrate and operate your weighing equipment with security
- Comply with quality and compliance standards in laboratory and manufacturing

www.mt.com/GWP

			/	ietura
1A/1A/1A/	mı	COM	/mn	ICTIILD

For more information

Mettler-Toledo GmbH

Im Langacher 44 8606 Greifensee, Switzerland www.mt.com/contact

Subject to technical changes.
© Mettler-Toledo GmbH 01/2021 30258547J en



30258547