# METTLER TOLEDO

## DGI117-Water

**Glass electrode with Sensor Chip**

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## 1. Start-up maintenance

- 1.1 The internal buffer solution (3) must cover the whole inside of the glass membrane (1). Remove air bubbles by gently shaking in vertical direction.
- 1.2 To equalize pressure remove rubber cap (8) or perforate its nipple (9).
- 1.3 Refill the bridge electrolyte (7) – (KCl 1 mol/L, PN 51343181)– to approx. 1 cm below the aperture (8).
- 1.4 To activate the glass membrane place the electrode in a tenfold diluted bridge electrolyte solution. Use a 0.5 mol/L KCl solution if no bridge electrolyte is at hand.

## 2. Plug & Play performance

The Sensor Chip which stores the sensor parameter is located in the screw cap of the sensor.

The sensor is recognized after connection to the sensor input of the titrator as well as after each restart of the system by reading the sensor parameter on the chip. A popup appears on the titrator terminal when the sensor was successfully recognized. This popup allows you to directly enter the sensor setup. You find further information concerning the functionality of the Plug & Play sensor.

## 3. Adjustment and calibration

The parameters zero-point and slope are measured by adjustment. A new electrode should have parameter values according to 3.1–3.2. Instructions for the adjustment method are given in the titrator manual.

- 3.1 The zero point of the electrode (pH value at  $E = 0$  mV) is approx. pH 7. The electrode potential, measured in a buffer solution of pH 7, must be within  $-30$  mV and  $+30$  mV.
- 3.2 The electrode slope (potential change per pH unit) at the pH range 4–9 must be greater than  $-55$  mV/pH (at  $25$  °C).
- 3.3 You should test the response time of the electrode. After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 3.4 If calibrating a used electrode the zero point should be within pH 6 and 8, the slope not less than  $-52$  mV/pH, and the potential change after 60 s should not exceed 3 mV/30 s.
- 3.5 The correct operation of the Pt1000 temperature sensor is verified by measuring the temperature of ice water, which shall yield a value of  $0.0$  °C  $\pm 1$  °C. The temperature measurement is performed with the manual operation "Measure" in the titrator, which is described in the operating instructions of the titrator.

#### 4. Storage and lifetime

- 4.1 Store the electrode in the bridge electrolyte – the sleeve junction (4) has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 4.2 Never let the glass membrane dry out, otherwise it has to be reactivated and the KCl crystals in /on the sleeve junction have to be dissolved.
- 4.3 The lifetime of the electrode is 6 months to 3 years, depending on its use and treatment.

#### 5. Maintenance instructions

- 5.1 The measuring range extends from pH 1 to 11, the tolerable temperature range from 0 to 100 °C (32 to 212 °F).
- 5.2 The sample solution should never enter through the sleeve junction into the bridge electrolyte. Therefore the level of the bridge electrolyte must always be higher than the sample solution level.
- 5.3 Just wipe the electrode after rinsing, never rub it dry.
- 5.4 Do not let the cable connector plug corrode.
- 5.5 A dried out glass membrane can be reactivated by placing it for several hours in 0.1 mol/L HCl.

#### 6. Cleaning procedures

- 6.1 Contamination of the bridge electrolyte: Drain the bridge electrolyte through aperture (8) and proceed according to 1.3 and 1.4.
- 6.2 Blocked sleeve junction by AgCl: place it overnight in concentrated ammonia, rinse and condition it for 1 hour in a buffer solution pH 4.
- 6.3 Blocked sleeve junction by sulfide precipitate: place it in a solution of thiourea (8% thiourea/1 mol/L HCl, e.g. METTLER TOLEDO PN 51350102) until the dark color has vanished.
- 6.4 Blocked sleeve junction by protein: place it at least for 1 hour in pepsin solution (5% pepsin/0.1 mol/L HCl, e.g. METTLER TOLEDO PN 51350100).
- 6.5 Other contaminations: place it for a few minutes in an ultrasonic bath with deion. water, ethanol or a mixed acid solution (e.g. aqua regia).

#### 7. Regeneration

If the electrode does not meet the qualifications according to 3.4 anymore, try a corrosion of the glass membrane: place it 1–2 minutes in diluted HF solution (2% HF/5% HCl, e.g. METTLER TOLEDO PN 51350104), rinse thoroughly with deion. water and treat it as in 6.1.

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