# METTLER TOLEDO

## DMi140-SC/DMi144-SC

**Platinum ring electrode with Sensor Chip**

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**Plug & Play combined platinum ring electrode**  
with ceramic frit for redox titrations.

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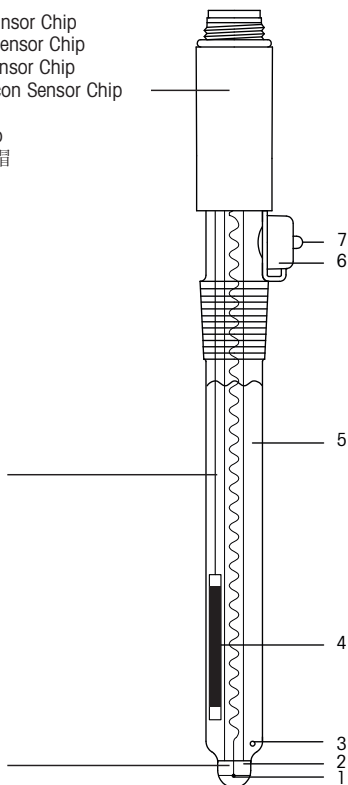
## 1. Start-up maintenance

- 1.1 To equalize pressure remove the rubber cap (6) or perforate its nipple (7).
- 1.2 Refill the reference electrolyte (5) – (e.g. METTLER TOLEDO No. 51340049) – to approx. 1 cm below the aperture (6).
- 1.3 Place the electrode in deion. water for 15 minutes to dissolve the KCl crystals in or on the frit (3).

## 2. Plug & Play performance

The Sensor Chip which stores the sensor parameter is located in the screw cap of the sensor.

The sensor is recognized after connection to the sensor input of the titrator as well as after each restart of the system by reading the sensor parameter on the chip. A popup appears on the titrator terminal when the sensor was successfully recognized. This popup allows you to directly enter the sensor setup. You find further information concerning the functionality of the Plug & Play sensor.

## 3. Testing

The electrode can be tested with a redox buffer solution, i.e. its potential and response time are measured.

- 3.1 The electrode potential should correspond to the values indicated on the buffer solution (e.g. METTLER TOLEDO No. 51340065 or 5134006).
- 3.2 Response time of the electrode: After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 3.3 Testing a used electrode, a comparison between titration curves is suitable: perform a titration with sodium thiosulfate (0.1 mol/L) and iodine (0.1 mol/L) with the new electrode. The titration goes fast with a big steep potential change at the equivalence point. A used electrode performing the same titration should present both these parameters.

## 4. Storage and lifetime

- 4.1 Store the electrode in the reference electrolyte – the frit has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 4.2 Never let the electrode go dry, otherwise the KCl crystals in or on the frit have to be dissolved.
- 4.3 The lifetime of the electrode is, depending on its proper use and treatment, almost unrestricted.

## 5. Maintenance instructions

- 5.1 The tolerable temperature range extends from 0 to 80 °C (32 – 176 °F).
- 5.2 In strong oxidizing solutions an oxygen adsorption takes place on the platinum surface of the electrode, in strong reducing solutions hydrogen adsorption. This can cause a change in the redox potential as well as an increased response time.
- 5.3 The sample solution should never enter into the reference electrode through the frit. Therefore the level of the reference electrolyte must be always higher than the sample solution level.
- 5.4 No reaction should take place between the sample solution and the reference electrolyte: sulfides, bromides, iodides and cyanides react with  $\text{Ag}^+$  and some cations ( $\text{Ag}^+$ ,  $\text{Hg}^{2+}$ ,  $\text{Au}^{3+}$ ,  $\text{Pb}^{2+}$ ) react with  $\text{Cl}^-$  of the reference electrolyte and cause a blockage of the frit.
- 5.5 Solutions containing protein quickly block the frit, due to protein precipitates.
- 5.6 Do not let the cable connector plug corrode.

## 6. Cleaning procedures

- 6.1 Contamination of the electrolyte: drain the electrolyte solution and refill with new electrolyte. Place the electrode in deion. water for several hours.
- 6.2 Frit blocked by  $\text{AgCl}$ : place it overnight in concentrated ammonia, rinse and condition it for 1 hour in a buffer solution pH 4.
- 6.3 Frit blocked by sulfide precipitate: place it in a solution of thiourea (8% thiourea /1 mol/ L HCl, e. g. METTLER TOLEDO No. 51340070) until the dark color has vanished.
- 6.4 Frit blocked by protein: place it for at least 1 hour in pepsin solution (5% pepsin /0.1 mol/L HCl, e.g. METTLER TOLEDO No. 51340068).
- 6.4 Frit blocked by protein: place it for at least 1 hour in pepsin solution (5% pepsin/0.1 mol/L HCl, e.g. METTLER TOLEDO No. 51340068).
- 6.5 Contamination of the platinum surface: depending on the contamination several cleaning procedures can be carried out.
  - a) Place the electrode for a few minutes in an ultrasonic bath with deion. water or ethanol.
  - b) Place the electrode for one minute in chromium sulfuric acid or aqua regia and rinse it thoroughly with deion. water.
  - c) Place the electrode for 30 minutes in a pH 4 buffer solution (Potassium hydrogen phthalate) saturated with quinhydrone to remove the absorbed oxygen or hydrogen.
  - d) Polish the platinum surface with  $\text{Al}_2\text{O}_3$  powder.

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