Puncture Knife for electrode LoT 406-M6-DXK-S7/25
Einstichmesser zu Elektrode LoT 406-M6-DXK-S7/25
Couteau de perforation pour électrode LoT 406-M6-DXK-S7/25


Order no. 59 900 386

Mettler-Toledo GmbH, Process Analytics
 Im Hackacker 15, CH-8902 Urdorf, Switzerland
 Phone +41 44 729 62 11

www.mt.com/pro

Subject to technical changes
 © 08/2020 METTLER TOLEDO. All rights reserved
 Printed in Switzerland. 59 906 009 B

For addresses of METTLER TOLEDO Market Organizations please go to: www.mt.com/pro-MOs

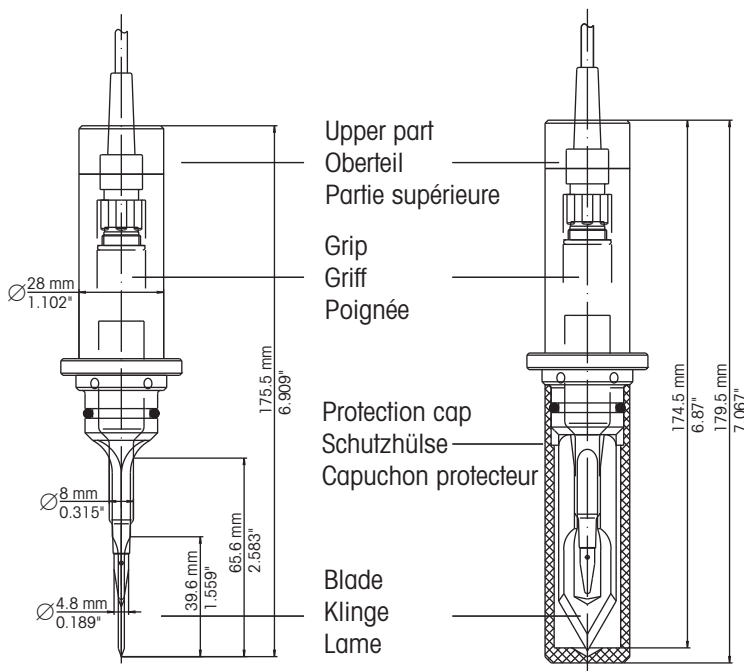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see up-side

2. Use

The puncture knife together with the electrode LoT 406-M6-DXK-S7/25 (Order no. 59 903 311), is used for direct pH measurement in meat, eliminating the need for pre-puncture.

3. Important notes

Store the knife in its protection cap!

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4. Instructions

4.1 Mounting

- a) Remove the watering cap from the electrode.
- b) Connect the cable to the electrode.
- c) Unscrew the upper part and the grip of the puncture knife.
- d) Slide the cable through the slit of the upper part and slide the grip over the electrode.
- e) Insert the electrode into the puncture knife (note position of the diaphragm aperture!).
- f) Screw the upper part and the grip onto. Choose the separate upper part provided when using the S7-BNC cable.
- g) Connect the cable to pH meter.

4.2 pH Measurement

Refer to the operating instructions for both the pH electrode and pH meter.

- Remove the protection cap before pH measurement.
- Calibrate (see pH electrode and pH meter instructions).
- Insert the puncture knife 3.5–6.5 cm into the medium and read the pH value.

4.3 Cleaning

The individual parts of the puncture knife can be cleaned with a commercially available detergent and a brush, followed by rinsing with ethanol.

4.4 Storage of the pH electrode in the puncture knife

- Place a little FRISCOLYT-B (Order no. 51 350 076) or 3 mol/L KCl (Order no. 51 350 072) in the protection cap of the knife.
- Replace the protection cap.

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