Operating Instructions

METTLER TOLEDO

Test Head QA000-W for the Quantos QB1 "MinWeigh" function

Introduction

The "MinWeigh" function defines a lower limit for the sample weight (in reference to the measurement uncertainty of the balance) in order to meet the admissible tolerance for the measured values, as specified by the work standard used (example: when working according to GMP the admissible tolerance is 1%, 2s, while the more restrictive USP standard allows for 0.1%, 3s). The "MinWeigh" function must be programmed and enabled and by a service engineer. In addition, the "MinWeigh" function must be activated in the menu in order to use the function (see Operating Instructions of the dosing system).

The QA000-W "MinWeigh" Test Head may be used by customers and METTLER-TOLEDO service engineers:

- Customers may use the test head to verify whether their dosing system still meets the requirements in terms of the minimum sample size (described below).
- METTLER-TOLEDO service engineers may use the test head to determine the minimum sample size on site, and to
 program the "MinWeigh" function accordingly (the procedure is described in the Service Manual).

Hardware requirements

- Quantos Powder Dosing System QB1 with terminal version V2.0 and later
- QA000-W Test Head
- Protocol printer RS-P25/QB1

The above configuration requires an Ethernet interface connected to an Ethernet-to-Serial converter (in most cases these items are part of the standard scope of delivery).

Preparation work

- Remove the test head from the storage container (same plastic container as used with standard dosing heads).
- Undo the knurled screw (C) at the top of the black barrel. The barrel holds the tare weight support (D) and 6 tare weights (A) of approximately 10g each. Take out the required number of tare weights making sure their total weight does not exceed the preset tare value of the "MinWeigh" function (see Operating Instructions of the dosing system, Chapter 6.13). When calculating the weight please also consider the tare weight support (D) which weighs in at approximately 5g. The overall tare (tare weights + support) must be slightly less than the programmed tare value!

Example: If the programmed tare value is 30g, the test should be carried out using a tare of approximately 25g which equals two tare weights (2 x 10g) plus the support (5g).

 Make sure the dosing system is operational and the "MinWeigh" function is active (see Operating Instructions of the dosing system).

Subject to technical changes and to changes in the accessories supplied with the instruments.

© Mettler-Toledo AG 2009 11781316 Printed in Switzerland 0904/2.32 Mettler-Toledo AG, Laboratory & Weighing Technologies, CH-8606 Greifensee, Switzerland

Test procedure

- Install the test head in the dosing unit. The dosing system then reads the head data.
- Make sure the appropriate "MinWeigh" method is activated (see Operating Instructions of the dosing system).

- Place the support (D) holding the tare weights (A) in the center of the weighing pan.
- Press the «Start» function key to launch the test procedure. You are now asked to adjust the test appliance.
- Turn the dosing head height adjustment handle to lower the dosing head until the tapered pin with the test weight (B) fully enters the pinhole of the tare weight support (D). This centers the tare weight support.
- Now turn the height adjustment handle back until the test weight is approximately 1 mm above the pinhole of the tare weight support (see figure).
- Using the arrow keys of the display you may move the test weight up and down in order to verify correct positioning of the tare weight support: in both end positions the display must show a stable weight value, otherwise you need to re-adjust the entire test appliance as described above.
- Press «OK» to confirm correct positioning of the test appliance.
- You are now asked to enter your User ID. Confirm your entry with «OK». This will start the test measurements.

Test measurements are fully automatic and the results are continuously sent to the protocol printer (in addition, the results are also transmitted to the host computer in XML format). After all measurements have been carried out, the final result will be printed:

- "PASSED": The test was successful and the programmed minimum sample size is still valid.
- "NOT PASSED": The test failed. Repeat the entire test procedure (including the adjustment of the test appliance). If the test fails several times it must be assumed that the programmed minimum sample size is no longer valid. In this case please call for a Mettler-Toledo service engineer.

Note:

The "LabXdirect quantos" software version 1.0 cannot process the XML file sent to the host computer. This feature will be implemented in a later version.

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