InSUS 300(i) Series

INGOLD
Leading Process Analytics

English	Installation Manual
Čeština	Montážní příručka
Dansk	Installationsvejledning
Deutsch	Installationshandbuch
Español	Manual de instalación
Suomi	Asennusopas
Français	Manuel d'installation
Italiano	Manuale di installazione
日本語	インストールマニュアル
한국어	설치 매뉴얼
Nederlands	Installatiehandleiding
Polski	Instrukcja instalacji
Português	Manual de Instalação
Svenska	Installationsmanual
ภาษาไทย	คู่มือการติดตั้ง
简体中文	安装手册

InSUS 30X(i)

InSUS 31X(i)

METTLER TOLEDO

Installation Manual	English
Montážní příručka	Čeština
Installationsvejledning	Dansk
Installationshandbuch	Deutsch
Manual de instalación	Español
Asennusopas	Suomi
Manuel d'installation	Français
Manuale di installazione	Italiano
インストールマニュアル	日本語
설치 매뉴얼	한국어
Installatiehandleiding	Nederlands
Instrukcja instalacji	Polski
Manual de Instalação	Português

Installationsmanual

Svenska

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Table of Contents

1	Introduction	3
2	Safety Instructions	3
3	Product Description	3
4	Installation	4
4.1	Unpacking.....	4
4.2	Installation into an Eldon James bag port	4
5	Storage	6
6	Environmental Protection	6
7	Warranty	6

1 Introduction

Thank you for buying this METTLER TOLEDO product. InSUS™ pH sensors are gamma and X-ray irradiation-sterilizable, pre-calibrated single-use pH sensors with an integrated Pt 1000 temperature probe. The sensors have been designed especially for deployment in single-use bags used in biotechnical processes. The sensors offer a design that allows safe and easy installation in a standard Eldon James port disc with a 1" barb. This port disc is welded into the bag film prior to sensor installation.

For installation into any port not described in this manual, contact the port manufacturer for instructions.

2 Safety Instructions

- The sensor contains glass components and should be handled carefully.
- Make sure that installation and operation is carried out only by trained personnel who have read and understood the instructions.
- A defective product must neither be installed nor put into service.
- No modifications to the product and the accessories are allowed. The manufacturer/supplier accepts no responsibility for damage caused by unauthorized modifications. The risk is borne entirely by the user.

3 Product Description

1. VarioPin (VP) connector (without protective cap)	4. Sensor shaft (HDPE)
2. Sensor label (serial number, factory calibration data)	5. O-rings (EPDM)
3. Retainer ring	6. Glass elements (reference electrode, pH electrode, and temperature probe)

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Sensor Quality Certificate showing factory end control data.

4 Installation

4.1 Unpacking

On unpacking, check the sensor for possible mechanical damage of the glass elements, the O-rings and the Vario- Pin (VP) connector. Report any damage to your METTLER TOLEDO supplier.

4.2 Installation into an Eldon James bag port

CAUTION

Do not clean the sensor glass elements with water, solvents or any other cleaning tools!

CAUTION

Removal or partial destruction of the protective Kollidon® (PVP) coating on the glass reference electrode of an InSUS 3XX(i) sensors with a shelf life longer than 12 months must be prevented during sensor handling!

NOTICE

1" Eldon James ports are available in different polymer materials. For a leak proof sensor installation METTLER TOLEDO offers InSUS sensors with different O-ring configurations. If in doubt please contact your METTLER TOLEDO supplier.

Recommended port / sensor combinations are:

1" Eldon James port material	Compatible InSUS pH sensor types
Engage™ 8402	InSUS 30X(i), InSUS 31X(i)
Dowlex™ 2517	InSUS 30X(i)/S, InSUS 31x(i)/S

- If applicable in your production environment use a mixture of 70 % isopropyl alcohol and 30 % deionized water for carefully lubricating the sensor O-rings prior to the insertion into the Eldon James port.
- 1 Remove the retainer ring from the sensor.
 - 2 Push the sensor all the way into the port until the sensor flange is flush with the port rim.
 - ➔ This is ideally done with a slight twisting motion to facilitate the insertion.

**NOTICE**

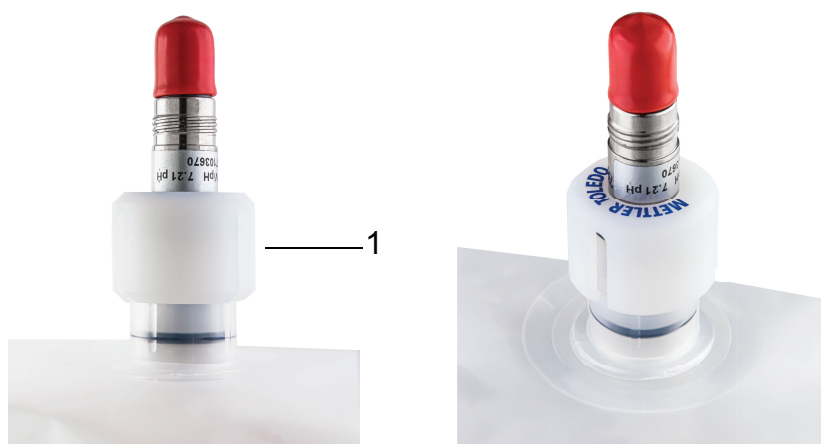
Make sure that the O-rings are still located completely in their grooves after this insertion procedure. If this is not the case, pull the sensor out of the port, reposition the O-rings and repeat the insertion procedure.

Correct sensor and O-ring positions after sensor insertion into an Eldon James port.

1. Shoulder	3. Port rim
2. Sensor flange	4. O-rings

- 3 Push the retainer ring over the sensor flange and port shoulder until it clicks into place.

Correct position of the retainer ring on sensor and bag port

1. Retainer ring

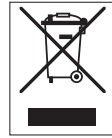
5 Storage

Storage conditions: Dry, 5 to 25 °C (41 to 77 °F), ambient pressure

6 Environmental Protection

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties, the content of this regulation must also be related.

7 Warranty

In the event of manufacturing faults, 12 months warranty from date of production is granted.

- InSUS is a trademark of the METTLER TOLEDO Group.
- Engage and Dowlex are trademarks of the Dow Chemical Company.
- Kollidon is trademark of BASF

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© 02/2026 METTLER TOLEDO. All rights reserved.
Mettler-Toledo GmbH