

for Analytical Balances XPE, XSE



Choose a place for the Smart-Tag

- Select a flat area (if only round areas are available, choose a bigger radius)
- Select a place without metal behind the sticking area
- Select a place outside the gripping area
- Select a place that is easy to hold to the reader.

Sticking the Smart-Tag to the pipette

1. Clean the sticking area with a mild solvent (e.g. Isopropanol, Metanol).
2. Wait until the area is completely dry.
3. Press the Smart-Tag to the area.
4. Leave it untouched for one day.

www.mt.com/excellence

For more information

Mettler-Toledo AG

Laboratory & Weighing Technologies
CH-8606 Greifensee, Switzerland

Subject to technical changes
© Mettler-Toledo AG 10/2013
30104632A en

