
Analytical Balances and Comparators

XPR

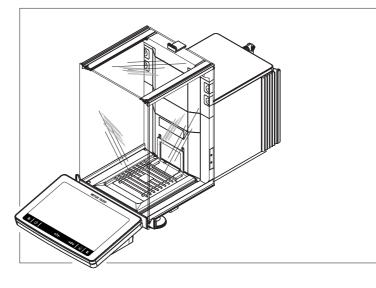






This User Manual provides brief instructions about the first steps to take with the instrument. This ensures safe and efficient handling. Personnel must have carefully read and understood this manual before performing any task.

For full information, always refer to the Reference Manual (RM).


www.mt.com/XPR-analytical-RM

Overview balance

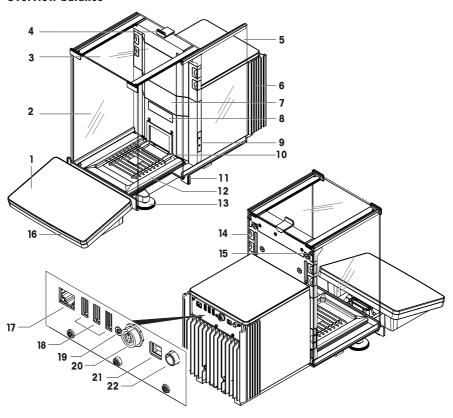

1	Terminal	12	Drip tray
2	Front panel draft shield	13	Leveling feet
3	Top door draft shield	14	Removable clips
4	Handle for top door	15	Side door release lever
5	Side door draft shield (right/left)	16	StatusLight
6	Cooling unit	17	Ethernet port
7	Slot for the installation of an internal module e.g., ionizer module	18	USB-A ports (to device)
8	Model label	19	Service seal
9	Optical sensor SmartSens	20	Socket for terminal connection cable
10	SmartGrid weighing pan	21	USB-B port (to host)
11	Door handle	22	Socket for AC/DC adapter

Overview terminal

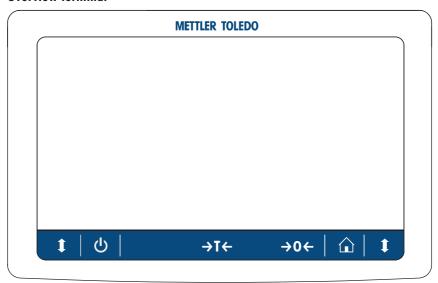

ψ	Standby	→T←	Tare
$\hat{\mathbf{u}}$	Home screen	→0 ←	Zero
‡	Open/close door		

Table of Contents

1	Introd	duction		3	
	1.1	Further do	ocuments and information	3	
	1.2	Acronyms	and abbreviations	3	
	1.3	Complian	ce information	4	
2 Safety Information					
	2.1		s of signal words and warning symbols	4	
	2.2		pecific safety information		
3	Desig	n and Fund	tions	5	
	3.1	Overview		5	
	3.2	User inter	face	5	
		3.2.1	Main sections at a glance		
		3.2.2	Main weighing screen	- 6	
4	Instal	lation and	Putting into Operation	7	
	4.1	Selecting	the location		
	4.2		g the balance		
	4.3	•	delivery		
	4.4		n		
		4.4.1	Attaching the terminal		
	4.5	4.4.2	Assembling the balance		
	4.5		to operation		
		4.5.1 4.5.2	Connecting the balance		
		4.5.2	Switching on the balance		
		4.5.4	Performing an internal adjustment		
		4.5.5	Entering / Exiting standby mode		
		4.5.6	Switching off the balance		
	4.6		g a simple weighing		
	1.0	4.6.1	Opening and closing the draft shield doors		
		4.6.2	Zeroing the balance		
		4.6.3	Taring the balance		
		4.6.4	Performing a weighing		
		4.6.5	Completing the weighing	15	
	4.7	Transport	ing, packing and storing	15	
		4.7.1	Transporting the balance over short distances	15	
		4.7.2	Transporting the balance over long distances	16	
		4.7.3	Packing and storing	16	
5	Main	tenance		16	
	5.1	Maintena	nce tasks	16	
	5.2	Cleaning		17	
		5.2.1	Disassembling for cleaning		
		5.2.2	Cleaning the balance		
		5.2.3	Putting into operation after cleaning	19	
6	Techr	nical Data		19	
	6.1	General d	ata	19	
7	Dispo	sal		20	

1 Introduction

Thank you for choosing a METTLER TOLEDO balance. The balance combines high performance with ease of use.

Disclaimer for comparators

In this document, the term "balance" is used to describe both balances and comparators.

Comparators are characterized by their higher resolution compared to balances. They are mainly used for differential weighing applications, such as the calibration of standard weights. Beside standard balance tests, comparators have also been tested with differential repeatability (ABA repeatability) during production.

FIII A

The software in this product is licensed under the METTLER TOLEDO End User License Agreement (EULA) for Software

When using this product you agree to the terms of the EULA.

www.mt.com/EULA

1.1 Further documents and information

This document is available in other languages online.

Product page:

www.mt.com/XPR-analytical

Instructions for cleaning a balance, "8 Steps to a Clean Balance":

www.mt.com/lab-cleaning-guide

Search for software:

www.mt.com/labweighing-software-download

Explanation

Search for documents:

www.mt.com/library

For further questions, please contact your authorized METTLER TOLEDO dealer or service representative.

www.mt.com/contact

Original term

1.2 Acronyms and abbreviations

AC	Alternating Current
ASTM	American Society for Testing and Materials
DC	Direct Current
EMC	Electromagnetic Compatibility
FCC	Federal Communications Commission
GWP	Good Weighing Practice
HID	Human Interaction Device
ID	Identification
LED	Light-Emitting Diode
LPS	Limited Power Source
MAC	Media Access Control
MT-SICS	METTLER TOLEDO Standard Interface Command Set
NA	Not Applicable
OIML	Organisation Internationale de Métrologie Légale
	(International Organization of Legal Metrology)
RAM	Random Access Memory
RFID	Radio-frequency identification
RM	Reference Manual

SELV Safety Extra Low Voltage
SOP Standard Operating Procedure
SQC Statistical Quality Control

UM User Manual

USB Universal Serial Bus

USP United States Pharmacopeia

1.3 Compliance information

National approval documents, e.g., the FCC Supplier Declaration of Conformity, are available online and/or included in the packaging.

www.mt.com/ComplianceSearch



For further information, consult the Reference Manual (RM).

www.mt.com/XPR-analytical-RM

2 Safety Information

Two documents named "User Manual" and "Reference Manual" are available for this instrument.

- The User Manual is printed and delivered with the instrument.
- The electronic Reference Manual contains a full description of the instrument and its use.
- Keep both documents for future reference.
- Include both documents if you transfer the instrument to other parties.

Only use the instrument according to the User Manual and the Reference Manual. If you do not use the instrument according to these documents or if the instrument is modified, the safety of the instrument may be impaired and Mettler-Toledo GmbH assumes no liability.

2.1 Definitions of signal words and warning symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

Signal words

DANGER A hazardous situation with high risk, resulting in death or severe injury if not avoided.

WARNING A hazardous situation with medium risk, possibly resulting in death or severe injury if

not avoided.

CAUTION A hazardous situation with low risk, resulting in minor or moderate injury if not

avoided.

NOTICE A hazardous situation with low risk, resulting in damage to the instrument, other

material damage, malfunctions and erroneous results, or loss of data.

Warning symbols



General hazard



Notice

2.2 Product-specific safety information

Intended use

This instrument is designed to be used by trained staff. The instrument is intended for weighing purposes.

Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties.

Mettler-Toledo GmbH assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. Mettler-Toledo GmbH assumes that the instrument owner provides the necessary protective gear.

Safety notes



⚠ WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to death or injury.

- Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.



NOTICE

Damage to the instrument or malfunction due to the use of unsuitable parts

- Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

A list of spare parts and accessories can be found in the Reference Manual.

3 Design and Functions



For further information, consult the Reference Manual (RM).

www.mt.com/XPR-analytical-RM


3.1 Overview

See the sections "Overview" (graphics and legend) at the very beginning of this manual.


3.2 User interface

3.2.1 Main sections at a glance

The main weighing screen (1) is the central navigation point where all the menus and settings can be found. The **Balance menu** (2), **Methods** (3) and **Results** (4) open when tapping the drawers along the main weighing screen.

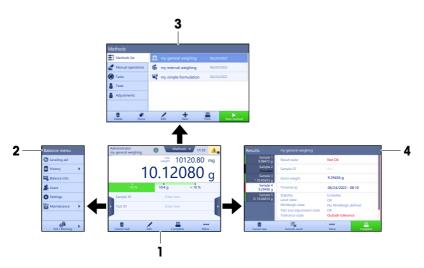

3.2.2 Main weighing screen

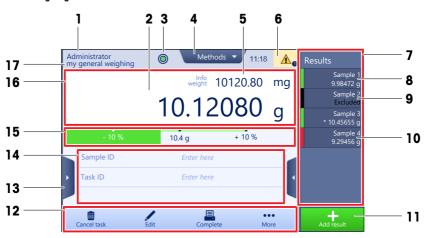

	Name	Description
1	User name	Shows the name of the current user.
2	Weighing value field	Shows the current weighing value.
3	Level indicator	Indicates if the balance is leveled (green) or not (red).
4	Methods menu	Accesses the user-defined list of methods, tests, and alignments.
5	Info weight	Shows the current weighing value in another unit.
6	Warning and error message area	Shows current warning and/or error messages.
7	Results list	Shows the weighing results saved for this task.

	Name	Description	
8	Sample status OK	Result status indicator green: indicates that the result fulfills a set of criteria. For example: The balance is in level. The internal adjustment was performed and ok. The weighing result is within the defined tolerance range (only if tolerance is defined).	
9	Sample status Excluded	Result status indicator black: indicates that the result was excluded from the Results list .	
10	Sample status Not OK	Result status indicator red: indicates that the result criteria are not fulfilled, e.g., "The weighing result was out of the defined tolerances".	
11	Button Add result	Adds the result to the Results list . Depending on the selected method, the button can have different functions.	
12	Action bar	Contains actions referring to the current task.	
13	Balance menu	Accesses the balance properties.	
14	Method information area	Contains information about the sample, method or task IDs.	
15	SmartTrac	Used as a weighing aid to define a target weight with upper and lower tolerances.	
16	Weighing value area	Shows the results of the current weighing process.	
17	Method name	Shows the name of the current method.	

4 Installation and Putting into Operation

4.1 Selecting the location

A balance is a sensitive precision instrument. The location where it is placed will have a profound effect on the accuracy of the weighing results.

Requirements of the location

Place indoors on stable Ensure sufficient spacing Level the instrument Provide adequate lighting table

Avoid direct sunlight Avoid vibrations Avoid strong drafts Avoid temperature fluctuations

Sufficient spacing for balances: > 15 cm all around the instrument Take into account the environmental conditions. See "Technical Data".


4.2 Unpacking the balance

Check the package, the packaging elements and the delivered components for damages. If any components are damaged, please contact your METTLER TOLEDO service representative.


i Note

Depending on the balance model, the components may look different. The procedure is always the same.


 Open the box and lift the package out using the lifting strap (1).

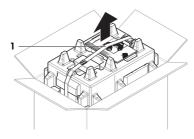


Open the lifting strap (1) and remove the User Manual (2).

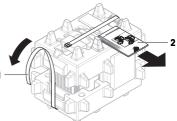


3 Remove the upper part of the package and remove the set with the AC adapter and power cable (3), the box containing several accessories (4), and the weighing pan (5).

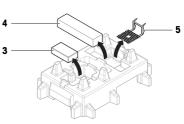


4 Carefully remove the terminal (6).

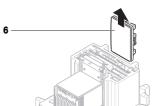

5 Carefully remove the package set with the draft shield doors and the display holder (7).

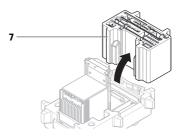

- 6 Carefully remove the platform (8) from the bottom packaging.
- 7 Remove the protective bag.
- 8 Store all parts of the packaging in a safe place for future
- → The platform is ready for assembling.

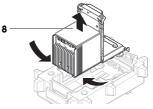

4.3 Scope of delivery

Balance

- Weighing unit
- Draft shield
- Drip tray and weighing pan

- Terminal with terminal holder and terminal connection cable
- AC/DC adapter with country-specific power cable
- MC Link Software (only comparators)

· Declaration of Conformity

Documentation

- User Manual
- Production certificate

Accessories

- ErgoClip basket 1,2
- SmartPrep², 2 pcs
- 1) not for XPR106DUH
- 2) not for XPR105DUHR and XPR106DUHR

Brush

4.4 Installation

4.4.1 Attaching the terminal




NOTICE


Damage to the cables due to careless handling

Do not kink or twist the cables.


1 Insert the slides of the display holder (1) into the front of the platform (2).

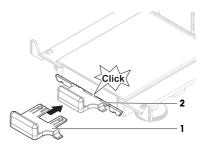


2 Connect the terminal cable (3) with the terminal (4). Consider the pin assignment.

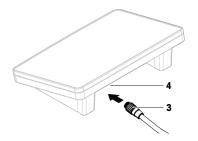


3 Place the terminal (5) onto the terminal holder (6).

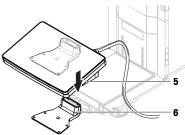

- 4 Carefully tilt the balance to its side.
- 5 Lead the cable (7) through the cable channel (8).
- 6 Carefully put the balance back on its feet.

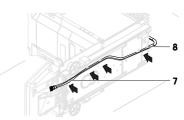

- 7 Insert the plug of the terminal cable (9) into the socket of the balance (10).
- → The terminal is ready.

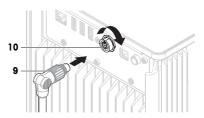

4.4.2 Assembling the balance



CAUTION

Injury due to sharp objects or broken glass


Instrument components, e.g., glass, can break and lead to injuries.

- Always proceed with focus and care.


i Note

Depending on the balance model, the components may look different. The procedure is always the same.


- 1 Insert the drip tray (1).
- 2 Carefully mount the weighing pan (2).

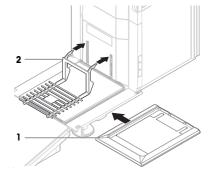


3 Place the side doors (3) into the grooves of the door slides (4) and tilt them up until they engage with the door lever (5). Consider the marks on the bottom frames (L = left / R = right).

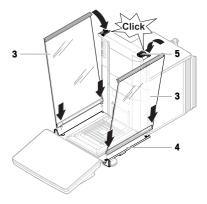

- 4 Insert the front panel (6) into the grooves (7) and tilt it up until it engages.
- 5 Open the side doors.

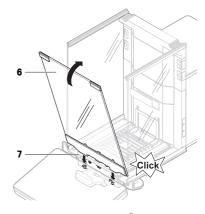

- 6 Fit the top door (8) along the top frame of the side doors and into the rails of the back wall (9).
- 7 Push the top door (8) towards the front.
- 8 Close the side doors.
- The balance is assembled and ready to be put into operation.

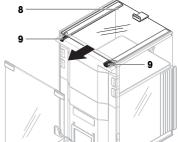

4.5 Putting into operation

4.5.1 Connecting the balance




↑ WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to death or injury.

- Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.
- 1 Install the cables in such a way that they cannot be damaged or interfere with operation.
- 2 Insert the plug of the AC/DC adapter (1) into the power socket of the instrument.
- 3 Secure the plug by firmly tightening the knurled nut.
- 4 Insert the plug of the power cable into a grounded power outlet that is easily accessible.
- ⇒ The balance automatically switches on.
- The draft shield opens and closes for initialization.

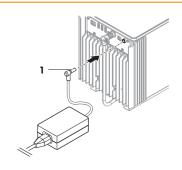

i Note

Do not connect the instrument to a power outlet controlled by a switch. After switching on the instrument, it must warm up before giving accurate results.

See also

General data ▶ Page 19

4.5.2 Switching on the balance

When connected to the power supply, the balance automatically switches on.

EULA (End User License Agreement)

When the balance is switched on the first time, the EULA (End User License Agreement) appears on the screen.

- 1 Read the conditions.
- 2 Tap I accept the terms in the license agreement, and confirm with \checkmark OK.
 - → The main weighing screen appears.

Acclimatization and warm up

Before the balance gives reliable results, it must:

- · acclimatize to the room temperature
- warm up by being connected to the power supply

The acclimatization time and warm-up time for balances and comparators are available in "General data".

Note

When the balance is exiting standby, it is ready immediately.

See also

- General data ▶ Page 19
- Entering / Exiting standby mode ▶ Page 14

4.5.3 Leveling the balance

Exact horizontal and stable positioning are essential for repeatable and accurate weighing results.

If the message Balance is out of level appears:

- 1 Tap ▶ Level the balance.
 - → The Leveling aid opens.
- 2 Turn both leveling feet (1) as instructed on the display until the dot is in the center of the level indicator.

The leveling aid can also be accessed through the **Balance** menu:

Navigation: ▶ Balance menu > ⊙ Leveling aid

4.5.4 Performing an internal adjustment

- The adjustment Strategy is set to Internal adjustment.
- 1 Open the Methods section, tap Adjustments, select the adjustment, and tap Start or -

from the main weighing screen, tap ... More and tap Start adjustment.

- → Internal adjustment is being executed.
- ⇒ When the adjustment has been completed, an overview of the adjustment results appears.
- 2 Tap **Print** if you want to print the results.
- 3 Tap **/ Finish adjustment**.
- → The balance is ready.

Navigation: ▼ Methods > • Adjustments

4.5.5 Entering / Exiting standby mode

- 1 To enter standby mode, hold **(b)**.
 - The display is dark. The balance is still switched on.
- 2 To exit standby mode, press **(b**).
 - → The display is turned on.

4.5.6 Switching off the balance

To completely switch off the balance, it must be disconnected from the power supply. By holding $oldsymbol{U}$, the balance goes only into standby mode.

i Note

When the balance was completely switched off for some time, it must warm up before it can be used.

See also


Switching on the balance ▶ Page 13

4.6 Performing a simple weighing

4.6.1 Opening and closing the draft shield doors

 Open the door manually with the door handle (1) or touch the key \$ on the terminal (2).

The doors can be configured to open and close in different ways.

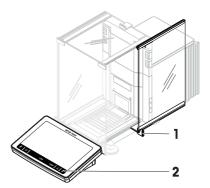


4.6.2 Zeroing the balance

- 1 Open the draft shield.
- 2 Clear the weighing pan.
- 3 Close the draft shield.
- 4 Press $\rightarrow 0 \leftarrow$ to zero the balance.
- ⇒ The balance is zeroed.

4.6.3 Taring the balance

If a sample vessel is used, the balance must be tared.

- 1 Open the draft shield.
- 2 Clear the weighing pan.
- 3 Close the draft shield.
- 4 Press $\rightarrow 0 \leftarrow$ to zero the balance.
- 5 Open the draft shield.
- 6 Place the sample vessel (1) on the weighing pan (2).
- 7 Close the draft shield.
- 8 Press \rightarrow **T** \leftarrow to tare the balance.
- → The balance is fared. The icon Net appears.

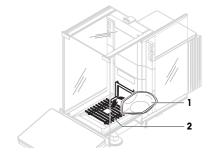

4.6.4 Performing a weighing

- 1 Open the draft shield.
- 2 Place the weighing object into the sample vessel.
- 3 Close the draft shield.
- 4 Tap + Add result if you want to report the weighing result.
- → The result is added to the **Results list**.

4.6.5 Completing the weighing

- 1 To save the **Results list**, tap **Complete**.
 - ⇒ The window Complete task opens.
- 2 Select an option to save or print the Results list.
 - → The respective dialog opens.
- 3 Follow the instructions from the wizard.
- 4 Tap Complete.
- The **Results list** is saved/printed and then cleared.


4.7 Transporting, packing and storing

4.7.1 Transporting the balance over short distances

- Disconnect the AC/DC adapter and unplug all interface cables.
- 2 Hold the platform with both hands and carry the balance in horizontal position to the target location. Consider the requirements of the location.

If you want put the balance into operation, proceed as follows:

- 1 Connect in reverse order.
- 2 Level the balance.
- 3 Perform an internal adjustment.

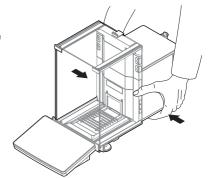

See also

- Selecting the location ▶ Page 7
- Switching on the balance ▶ Page 13

4.7.2 Transporting the balance over long distances

METTLER TOLEDO recommends using the original packaging for transportation or shipment of the balance or balance components over long distances. The elements of the original packaging are developed specifically for the balance and its components and ensure maximum protection during transportation.

See also

Unpacking the balance ▶ Page 8

4.7.3 Packing and storing

Packing the balance

Store all parts of packaging in a safe place. The elements of the original packaging are developed specifically for the balance and its components, and ensures maximum protection during transportation and storage.

Storing the balance

Only store the balance under the following conditions:

- · Indoor and in the original packaging
- · According to the environmental conditions, see "Technical Data"

i Note

When storing for longer than 6 months, the rechargeable battery may become empty (only date and time get lost).

See also

5 Maintenance

To guarantee the functionality of the balance and the accuracy of the weighing results, a number of maintenance actions must be performed by the user.



For further information, consult the Reference Manual (RM).

www.mt.com/XPR-analytical-RM

5.1 Maintenance tasks

Maintenance action	Recommended interval	Remarks
Performing an internal adjustment	DailyAfter cleaningAfter levelingAfter changing the location	see "Performing an internal adjustment"
Performing routine tests (eccentricity test, repeata- bility test, sensitivity test). METTLER TOLEDO recommends to at least perform a sensitivity test.	After cleaning After assembling the balance After a software update Depending on your internal regulations (SOP)	see "Tests" in the Reference Manual
Cleaning	After every use After changing the substance Depending on the degree of pollution Depending on your internal regulations (SOP)	see "Cleaning"

Maintenance action	Recommended interval	Remarks
Updating the software	 Depending on your internal regulations (SOP). After a new software release. 	see "Software update" in the Reference Manual

See also

- Performing an internal adjustment ▶ Page 13
- Cleaning ▶ Page 17

5.2 Cleaning

5.2.1 Disassembling for cleaning




CAUTION


Injury due to sharp objects or broken glass

Instrument components, e.g., glass, can break and lead to injuries.

- Always proceed with focus and care.
- 1 Open the top door (1) and pull it all the way back, outside of the rails of the side doors. Shortly before the top door drops out, you can feel a slight resistance. Just keep pulling a little bit tighter.

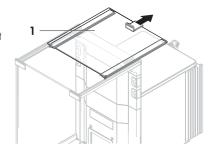


- 2 Hold the side doors (2) and push down the lever (3) to release them.
- 3 Carefully remove both side doors (2).

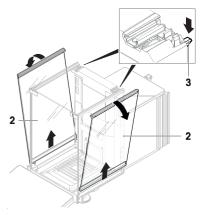

4 Tilt the front panel (4) to the front and remove it.



- 5 Carefully lift the weighing pan (5) to unhook it and pull it out
- 6 Remove the drip tray (6).
- 7 Store all removed components in a safe place.
- The balance is ready for cleaning.

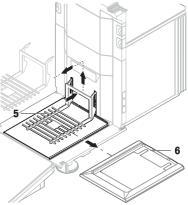

5.2.2 Cleaning the balance



NOTICE

Damage to the instrument due to inappropriate cleaning methods

If liquid enters the housing, it can damage the instrument. The surface of the instrument can be damaged by certain cleaning agents, solvents, or abrasives.

- 1 Do not spray or pour liquid on the instrument.
- 2 Only use the cleaning agents specified in the Reference Manual (RM) of the instrument or the guide "8 Steps to a Clean Balance".
- 3 Only use a lightly moistened, lint-free cloth or a tissue to clean the instrument.
- 4 Wipe off any spills immediately.



For further information on cleaning a balance, consult "8 Steps to a Clean Balance".

www.mt.com/lab-cleaning-guide

Cleaning around the balance

- Remove any dirt or dust around the balance and avoid further contaminations.

Cleaning the terminal

- Clean the terminal with a damp cloth or a tissue and a mild cleaning agent.

Cleaning the removable parts

 Clean the removed part with a damp cloth or a tissue and a mild cleaning agent or clean in a dishwasher up to 80 °C.

Cleaning the balance

- 1 Disconnect the balance from the AC/DC adapter.
- 2 Use a lint-free cloth moistened with a mild cleaning agent to clean the surface of the balance.
- 3 Remove powder or dust with a disposable tissue first.
- 4 Remove sticky substances with a damp lint-free cloth and a mild solvent, e.g., isopropanol or ethanol 70%.

5.2.3 Putting into operation after cleaning

- 1 Reassemble the balance.
- 2 Check that the draft shield doors (top, sides) open and close normally.
- 3 Check if the terminal is connected to the balance.
- 4 Reconnect the balance to the AC/DC adapter.
- 5 Check the level status, level the balance if necessary.
- 6 Respect the warm-up time specified in the "Technical Data".
- 7 Perform an internal adjustment.
- 8 Perform a routine test according to the internal regulations of your company. METTLER TOLEDO recommends performing a sensitivity test after cleaning the balance.
- 9 Press $\rightarrow 0 \leftarrow$ to zero the balance.
- → The balance is ready to be used.

See also

- Technical Data ▶ Page 19
- Performing an internal adjustment ▶ Page 13

6 Technical Data

6.1 General data

Power supply

AC/DC adapter (model no. Input: $100 - 240 \text{ V AC} \pm 10\%$, 50 - 60 Hz, 1.8 A

FSP060-DHAN3): Output: 12 V DC, 5 A, LPS, SELV

AC/DC adapter (model no. Input: $100 - 240 \text{ V AC} \pm 10\%$, 50 - 60 Hz, 1.5 A

FSP060-DIBAN2): Output: 12 V DC, 5 A, LPS, SELV
Cable for AC/DC adapter: 3-core, with country-specific plug

Balance power consumption: $12 \text{ V DC} \pm 10\%$, 2.25 A

Protection and standards

Overvoltage category: II
Degree of pollution: 2

Standards for safety and EMC: See Declaration of Conformity
Range of application: Use only indoors in dry locations

Environmental conditions

The limit values apply when the balance is used under the following environmental conditions:

Height above mean sea level: Up to 5000 m Ambient temperature: +10 - +30 °C

Temperature change, max.: 5 °C/h

Relative humidity: 30 – 70%, non-condensing

Acclimatization time: At least **8 hours** after placing the instrument in the same location

where it will be put into operation.

Warm-up time: At least 120 minutes after connecting the balance to the power

supply. When switched on from standby, the instrument is ready

for operation immediately.

The balance can be used under the following environmental conditions. However, the weighing performances of the balance may be outside the limit values:

Ambient temperature: $+5 \, ^{\circ}\text{C} - +40 \, ^{\circ}\text{C}$

Relative humidity: 20% to max. 80% at 31 °C, decreasing linearly to 50% at

40 °C, non-condensing

The balance can be disconnected and stored in its packaging under the following conditions:

Ambient temperature: $-25 - +70 \, ^{\circ}\text{C}$

Relative humidity: 10 - 90%, non-condensing

Environmental conditions for comparators

Comparators need to be used under the following environmental conditions to reach the specified performances:

Air speed, max.: 0.15 m/s

7 Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE), this equipment may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this equipment in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this equipment. Should this equipment be passed on to other parties, the content of this directive must also be passed on to the other party.



To protect your product's future:

METTLER TOLEDO Service assures the quality, measuring accuracy and preservation of value of this product for years to come.

Please request full details about our attractive terms of service.

www.mt.com/service

www.iiii.coiii/	excemence	-anaiyiicai

For more information

Mettler-Toledo GmbH

Im Langacher 44 8606 Greifensee, Switzerland www.mt.com/contact

Subject to technical changes.
© 11/2024 METTLER TOLEDO. All rights reserved.
30419872F en

