

Weight Verification System for Traceable Balance Testing



Proven Traceability

The built-in balance testing application records the weight-specific data and the test results. A detailed test report can be printed out with an external printer to provide full documentation of the performed balance test.



Enhanced Security

WeightLink™ will only allow the use of a valid test weight. Verification of the test weight prior to use means you can rest assured that your testing procedures are fully compliant.



Improved Productivity

Automatic data transfer of weight parameters into the balance firmware is much faster than manual data entry and removes the need to check and re-check handwritten entries. Printing out results is quick and easy.




Seamless and Secure Semi-Automated Weight Check

WeightLink™ is an innovative weight identification system that guarantees full traceability of your balance tests.

Each test weight has a unique identification number (UIN) on the base which 'links' it to its certificate. Calibration and certificate data is stored in a data matrix code (DMC). The dedicated DMC Scanner reads the two codes and sends the information to the balance. The balance firmware validates the certificate and weight data before allowing the balance test to begin.

With this highly efficient system, it is impossible to test using the wrong weight and the risk of errors from manual data entry is completely eliminated.

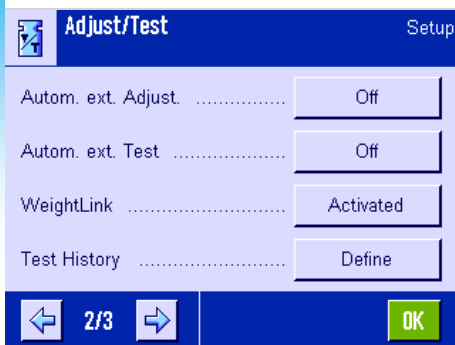

Easy Setup

WeightLink™ weights are available from 1 g to 5 kg and the data matrix scanner simply plugs into the back of your balance. Your local METTLER TOLEDO representative can advise you on compatible balance firmware versions.

The WeightLink™ System

WeightLink™ Weights

- Calibration certificate included
- Unique identification number on base
- Available individually or in a CarePac®

- Lifetime warranty and guaranteed positive tolerances*
- High grade stainless steel
- Supplied in an easy-clean plastic box

*Terms and Conditions can be found under www.mt.com/signature-line

WeightLink™ CarePacs®

Weight 1		Weight 2		
Nominal Weight / Class		Nominal Weight / Class		
200 g	F2	20 g	F1	30293475
200 g	F2	10 g	F1	30293476
100 g	F2	5 g	E2	30293477
50 g	F2	2 g	E2	30293478
2 g	E2	0.1 g	E2	30293479
5 g	E2	0.2 g	E2	30293480
20 g	F1	1 g	E2	30293481
500 g	F2	20 g	F1	30293482
1000 g	F2	50 g	F2	30293483
2000 g	F2	100 g	F2	30293484
2000 g	F2	200 g	F2	30293485
5000 g	F2	200 g	F2	30293486

Individual WeightLink™ Weights

Nominal weight	Class E2	Class F1
1 g	30293502	30293561
2 g	30293503	30293562
5 g	30293504	30293563
10 g	30293505	30293564
20 g	30293506	30293565
50 g	30293507	30293566
100 g	30293508	30293567
200 g	30293509	30293568
500 g	30293550	30293569
1 kg	30293551	30293570
2 kg	30293552	30293571
5 kg	30293553	30293572

WeightLink™ CarePac® Optional 3rd Weights

Nominal weight	Class E2
1 g	30293554
2 g	30293555
5 g	30293556
10 g	30293557
20 g	30293558
50 g	30293559
100 g	30293560



Our CarePacs® contain just two weights for testing at 5% and 100% of your balance capacity in accordance with USP <41> recommendations. An optional third weight can be selected for customer-specific testing requirements. To find the right CarePac® for your balance, please visit:

► www.mt.com/carepacs

Calibration eData



All necessary calibration data is stored in a data matrix code (eData) which is linked to the correct weight via the UIN. For convenient balance testing, the code is printed on the WeightLink™ Card or adhered to the weight box.


DMC Scanner



The scanner connects directly to your balance via the RS232 port. Image analysis technology is used to scan the UIN on the weight and the code with the calibration certificate eData.

WeightLink™ DMC Scanner	30268560
WeightLink™ DMC Scanner + RS232 option	30304696

Balance Firmware

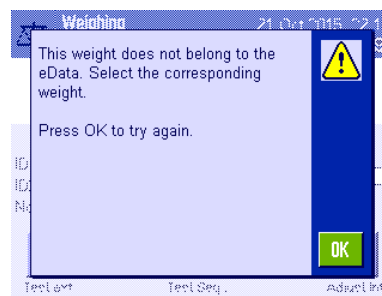

The balance firmware validates the certificate and weight data before allowing the balance test to begin. Please contact your local METTLER TOLEDO representative for more information on compatible balance firmware versions and their availability.



Mettler-Toledo GmbH

Laboratory Weighing
CH-8606 Greifensee, Switzerland
Phone +41 44 944 22 11
Fax +41 44 944 30 60

©11/2015 Mettler-Toledo GmbH
30303069
Global MarCom Switzerland, 1825 PH

www.mt.com/weightlink

For more information