



Project: BalanceLink V4.1.3	Product Documentation	
	Release Notes	Software Development Center (SWDC)

Release Notes METTLER TOLEDO BalanceLink		
Release: METTLER TOLEDO BalanceLink: 4.1.3	Release Date:	24. 07. 2019

Contents

1	Executable Signatures	2
2	Help File Signatures	2
3	Operating System Compatibility	2
4	System Hardware Recommendation	2
5	Modifications and solved issues.....	3
6	Notes	4

Project: BalanceLink V4.1.3	Product Documentation	
	Release Notes	Software Development Center (SWDC)

1 Executable Signatures

#	Filename	Version	Build	Date	Size [Bytes]
1	BAL32.exe	4.1.3	4.1.3.5	02.07.2019	2'002'344

2 Help File Signatures

Filename	Date		Filename	Date	
Bal32-E.chm	07.06.2019		Bal32-I.chm	07.06.2019	
Bal32-D.chm	07.06.2019		Bal32-S.chm	07.06.2019	
Bal32-F.chm	07.06.2019				

3 Operating System Compatibility


Software Description	Version	Service Packs	Remarks
Operating Systems (32/64 bit)	- Microsoft® Windows 7™	SP1	n/a
Operating Systems (32/64 bit)	- Microsoft® Windows 8.1™	None	n/a
Operating Systems (32/64 bit)	- Microsoft® Windows 10™	None	n/a

Notes:

Microsoft® (MS) Windows Operating Systems 95™ / 98™ / ME™ / NT™ / 2000™ / XP™ / Vista™ / 8.0™ are no longer supported/tested.


4 System Hardware Recommendation

Description	Min. System Requirements	Remarks
Processor	Any	
RAM	Any	
Hard Disk	20 MB available	
Screen Resolution	800 x 600	

Project: BalanceLink V4.1.3	Product Documentation	
	Release Notes	Software Development Center (SWDC)

5 Modifications and solved issues

Ticket	Description	Version
1	Supporting virtual COM ports (USB, IR)	V4.0.1
2	Password protection for configuration	V4.0.1
3	Supporting analytical instruments: SevenEasy S20 / S30; SevenGoPro	V4.0.1
4	Supporting moisture analyzers: HB43, HG63, HR83	V4.0.1
5	Supporting balances: AX, MX, UMX, XS, AB-S, PB-S, XP, SR, PL-S, Panda, BBA, BBK, IND	V4.0.1
6	Generic interface to connect not listed devices	V4.0.1
7	Language support for Spanish and Italian	V4.0.1
8	30 day trial version	V4.0.1
9	Now supporting MS Windows Vista™	V4.0.2
10	The Device Properties for Generic, S20, S30, SevenGoPro were not being saved properly to the BAL file.	V4.0.3
11	Trial version has no functional limitation anymore (formerly, there was e.g. no saving and no printing). Only the 30 day restriction remains.	V4.0.3
12	Support of the new big icons in explorer view under MS Windows Vista™ or Windows 7™ respectively.	V4.0.4
12	Support of MS Windows 8™. New CHM help files instead of HLP. New codec for movies on installation CD. Add-on of 1.2.2016: Compatibility tests were successfully done with MS Windows 8.1™ / Windows 10™ and MS Office 2010™ / MS Office 2013™/ MS Office 2016™	V4.1.0
13	Fix of issue while using Shift+Tab in Macro's	V4.1.1
14	Fix of Issue in 4.1.1 (only): If device property 'Do not modify (as is)' is used then there was a problem with an automatic 'Shift' add-on that led to strange highlighting and other effects within Windows Desktop.	V4.1.3
15	Support of USB driver for XPR/XSR balances.	V4.1.3
16	Updated 'Mettler-Toledo End User License Agreement (EULA) for Software' (and it is now in English for all languages).	V4.1.3
17	There is no section 'Moisture Analyzer' nor 'Instrument' available anymore. Those devices are supported by other MT software products. Please approach your sales representative.	V4.1.3
18	MS Windows Operating Systems 2000™ / XP™ / Vista™ / 8.0™ are no longer supported / tested.	V4.1.3
19	Compatibility tests were successfully done with MS Windows 10™ (version 1809) and MS Office 2019™.	V4.1.3

Project: BalanceLink V4.1.3	Product Documentation	
	Release Notes	Software Development Center (SWDC)

6 Notes

Installation:


Software must be installed with a supervisory-level login user name, such as "Administrator".

Executing the installation file a second time and confirming the 'Confirm Uninstall' dialog leads to the complete removal of all files installed except the configuration files.

Registration:

Registration must be done with a supervisory-level login user name, such as "Administrator".

Depending on the chosen directory it is sometimes necessary to run BAL32.exe once using right-click and 'Run as administrator' to make a successful registration:

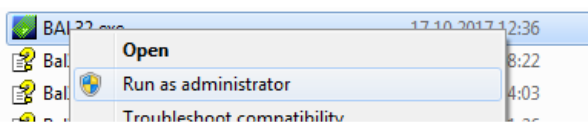

Add-On to Macro (View / Transmitter / Properties / Macro 1, 2, 3) - Supported Navigation Commands:

1. Cursor navigation combined commands (they are considered pressed together):

Ctrl+Up	
Ctrl+Left	
Ctrl+Home	(Cell A1)
Ctrl+End	(Furthest used cell)
Shift+Tab	

2. Cursor navigation commands:

Page Up
 Page Down
 Right
 Left
 Up
 Down
 Tab
 Home
 End
 Backspace
 Enter