USB Application: Basic, Count, Count+, Smart+

The 4-Series USB port is a USB slave intended for connection to a PC. It is not intended to connect to printers, scanners, etc. because the port is Slave, not Master.

In this demo we will connect a 4-Series Basic (BBA422) to a PC over USB and communicate bi-directional in dialog mode.

Setup the PC application.

We will demonstrate the application Hyper Terminal.

We determine the USB port assignment in our PC, here we use Windows XP.

Open Control Panel on the PC.

Click on "System"



On the System Properties tab	, click "Device Manager"
------------------------------	--------------------------

System Re	n Restore Automatic Updates Rem		Remote
General	Computer Name	Hardware	Advanced
	100		
Add Hardwa	re Wizard		
S 14	ie Add Hardware Wizard	i helps you install hards	ware.
~			
		Add Hardwar	e Wizard
NORMAN			
Hevice Man	ader		
Device Man	ager va Daviaa Managar lista	- 11 Maa la suduusus alsu üs	aa inatallad
Device Man	ager 1e Device Manager lists : 1 your computer. Use the	all the hardware device Device	es installed
Device Man Tł or pr	ager ne Device Manager lists 1 your computer. Use the operties of any device.	all the hardware device Device	es installed
Device Man Tł or pr	ager ne Device Manager lists i your computer. Use the operties of any device.	all the hardware device Device	es installed
Device Man Tł or pr	ager ne Device Manager lists i your computer. Use the operties of any device. Driver Signing	all the hardware device Device Device Ma	es installed the anager
Device Man Tł or pr	ager ne Device Manager lists i your computer. Use the operties of any device. Driver Signing	all the hardware device Device Device Ma	anager
Hardware Pr	ager he Device Manager lists hyour computer. Use the operties of any device. Driver Signing ofiles	all the hardware device Device Device Ma	anager
Hardware Pr	ager ne Device Manager lists your computer. Use the operties of any device. Driver Signing ofiles ardware profiles provide a ferent bardware configur	all the hardware device Device Ma Device Ma Device Ma	anager and store
Hardware Pr	ager ne Device Manager lists your computer. Use the operties of any device. Driver Signing ofiles ardware profiles provide a ferent hardware configur	all the hardware device Device Ma Device Ma a way for you to set up rations.	anager
Hardware Pr	ager he Device Manager lists your computer. Use the operties of any device. Driver Signing ofiles ardware profiles provide a ferent hardware configur	all the hardware device Device Ma Device Ma a way for you to set up rations.	anager
Hardware Pr	ager he Device Manager lists your computer. Use the operties of any device. Driver Signing ofiles ardware profiles provide a ferent hardware configur	all the hardware device Device Ma Device Ma a way for you to set up rations. Hardware	anager and store
Hardware Pr	ager ne Device Manager lists nyour computer. Use the operties of any device. Driver Signing ofiles ardware profiles provide a ferent hardware configur	all the hardware device Device Ma Device Ma a way for you to set up rations. Hardware	anager and store

On the Device Manager screen open "Ports (Com & LPT) to find where the USB port is configured (from the USB installation).



Here we find the USB (METTLER TOLEDO Serial Port) configured on COM 4 of our PC. With this information we can close all control panel windows and proceed with the Hyper Terminal application demo.

This demo is not on how to use Hyper Terminal program and a familiarity of it is assumed.

USB Properties		? 🗙
Connect To Set	tings	
See USB	Change Icon	
Country/region:	United States (1)	
Enter the area c	ode without the long-distance prefix.	
Area code:	1	
Phone number:		
Connect using:	COM4 Xircom Cardbus Ethernet 100 + Moder COM3 COM1	
Use country/	COM4 TCP/IP (Winsock) sy	
	ОК С	ancel

On the Hyper Term Properties Tab, select Connect to Com 4

On the Hyper Terminal "Connect To" tab, click the Configure button and set parameters to 9600 Baud, 8 data bits, no parity.

JSB Properties	? 🗙	
USB	COM4 Properties Port Settings	? 🗙
Country/region: United States (1) Enter the area code without the long- Area code: 1	Bits per second: Data bits:	9600
Phone number: Connect using: COM4 Configure	Parity: Stop bits:	None 💌
Use country/region code and area	Flow control:	Hardware 💌 Restore Defaults
		K Cancel Apply

Switch to the "Settings" tab and click the "ASCII Setup" button.

USB Properties	? 🛛
Connect To Settings	
 Function, arrow, and ct Terminal keys 	rl keys act as O Windows keys
Backspace key sends	O Ctrl+H, Space, Ctrl+H
Emulation:	
Auto detect	Y Terminal Setup
Telnet terminal ID:	ANSI
Backscroll buffer lines: 5	500 🗢
Play sound when con	necting or disconnecting
Input Translation	ASCII Setup
	OK Cancel

On the ASCII Setup page check mark the "Send line ends with line feeds". This is required for bidirectional communication with 4-Series.

ASCII Setup 🛛 💽 🔀
Sending Service in the ends with line feeds
Line delay: 0 milliseconds.
Character delay: 0 milliseconds.
ASCII Receiving Append line feeds to incoming line ends Force incoming data to 7-bit ASCII Wrap lines that exceed terminal width
OK Cancel

4 Series Scale Setup

In the setup of the 4-Series we will set COM 2 (because this is the USB port) to dialog mode and set the communication parameters.

Go into setup.

- Hold down the PRINT key until "CODE" is displayed
- Key in the Supervisor Password, then PRINT key.
- The top level menu will show SCALE, rotate through the menu with the "no" key, TARE until COMMUNICATIONS (CONNUNI) is displayed.
- Touch the "yes" key, PRINT, to enter the communications menu.
- COM 1 (CON I) is displayed. Rotate to COM 2 with the TARE key.
- Touch the "yes" key, PRINT, to enter the COM 2 menu.
- The mode (NOdE) prompt is displayed, touch the "yes" key, PRINT, to enter the mode menu.
- Rotate through the mode menu with the "no", TARE key until dialog is displayed (dIALOG), touch the "yes" key, PRINT to select dialog.

- The menu now advances to parameters (PArANEt). Touch the "yes" key PRINT.
- Set baud to 9600
- Set parity (Parity) to 8 none (8 nonE)
- Set handshake (HSHAkE) to NO or XONXOFF.
- Touch the ON/OFF key to short cut setup and exit.
- At the End and SAVE prompts touch the "yes" key, PRINT.

Run the demo using Hyper Terminal

Click the "Call" icon to start the session.

🌯 USB - HyperTermin	nal	
File Edit <u>View</u> Call Tr	ansfer Help	
0 🖻 🍘 🕐 🗈	ð 🕈	
Call		~
		~
<		>
Disconnected	Auto detect	9600 8

Now we can demonstrate communications by sending SICS commands and receiving the response from the 4-Series scale in the Hyper Terminal screen. Here are some sample SICS commands with the scale responses captured on the Hyper Terminal screen.

SI [Send weight immediate] I4 [Serial Number]

I2 [Scale model and capacity]



End of demo.