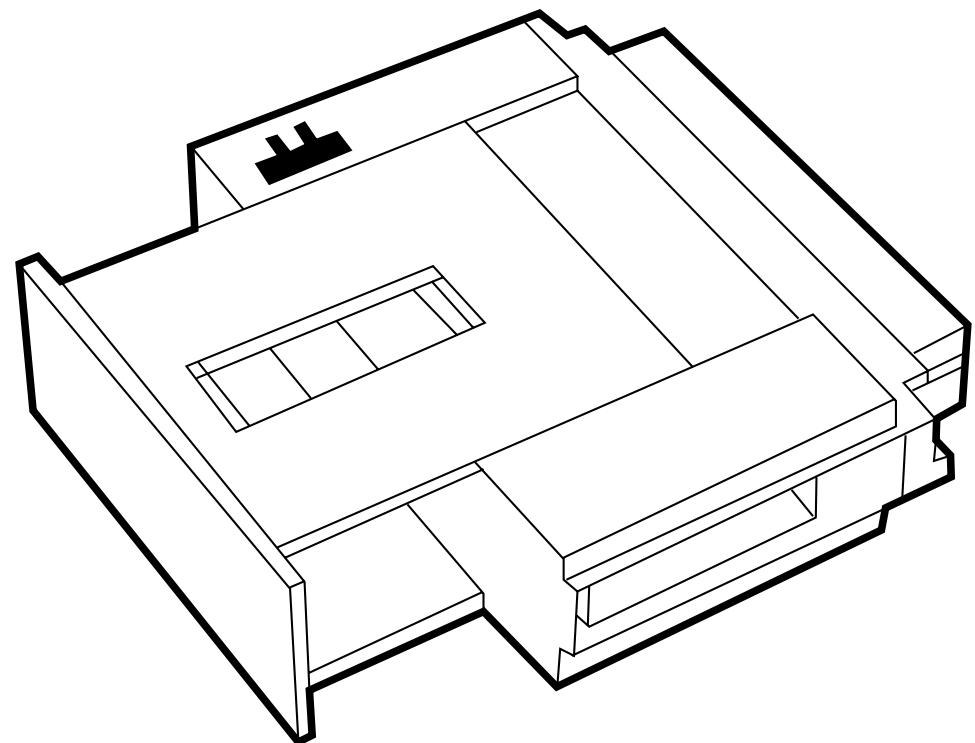
Operating Instructions
Bedienungsanleitung
Mode d'emploi
Istruzioni d'uso
Instrucciones de manejo

METTLER TOLEDO
CountEasy-M

METTLER TOLEDO

English Page 2

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
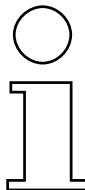
General

CountEasy-M is a program cassette for SM scales, high-capacity PM scales or for AM/PM balances up to 6.1 kg with the GM303 Control Unit. It allows you to use your balance or scale for convenient counting with a variable reference piece number.

The **insertion of the program cassette** is described in the operating instructions of the balance or scale in the section “Configuration”. When CountEasy-M is inserted, the application “Counting” can be activated with the **F1** key after the balance or scale has been switched on.

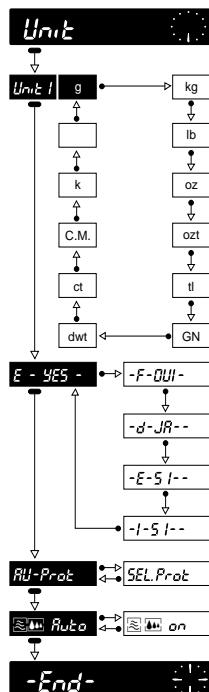Configuration register with CountEasy-M

The configuration register with CountEasy-M is distinguished from the configuration register of the balance or scale with the standard program cassette only in the –Unit– sector. With CountEasy-M the following can be selected in the –Unit– sector:

Language –E-YES– English (default setting)

–F-OUI– French

–d-JA– German

–E-SI– Spanish

–I-SI– Italian

Printout Au-Prot automatic printout (default setting)

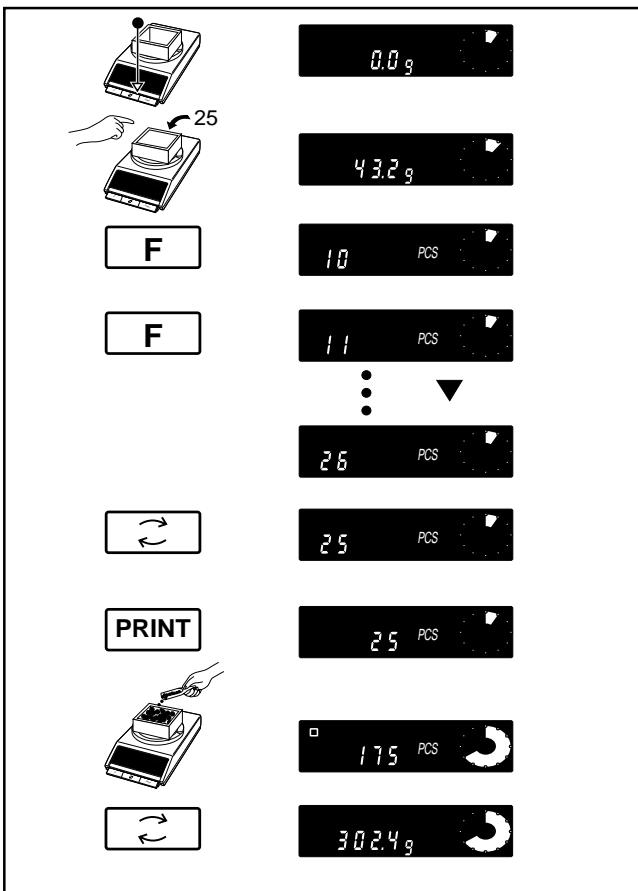
SEL.Prot printout only with the **PRINT** key

All further setting options are identical to those with the standard cassette, see operating instructions of the balance or scale.

Notes

- If the configuration register can not be selected, check whether the jumper of the program cassette is in the position “unsecured” (see operating instructions of the balance or scale, section “Configuration”).
- When the balance or scale is reset to the default configuration (Reset), the language is not reset.

Operation

Load empty or full container, tare.

Add reference piece number (25 Pcs) to the empty container or remove from the full container.

Recall of the set reference piece number: press key **F**.

This simultaneously switches from the weighing mode to the application "Counting".

Set reference piece number

- increment by 1: press key **F** briefly.
- increment continuously: press and hold key **F**.
- decrement by 1: press key **F** briefly.
- decrement continuously: press and hold key **F**.

The reference piece number is automatically stored if no key has been pressed for approx. 3 seconds or if during this time the **PRINT** key is pressed. Once the reference piece number has been stored, the calculated piece weight can be printed out with **PRINT**.

Add or remove (negative piece number) the number to be determined.


Switching between piece number and weight: press key **F**.

The piece number can be printed out with **PRINT**.

What if ...

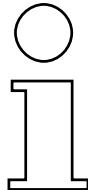- | |
|------------|
| Add |
|------------|
- Reference weight too low,
 - increase reference piece number (minimum piece weight = 1/4 digit).

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© Mettler-Toledo GmbH 2000 704712 Printed in Switzerland 0008/2.45
Mettler-Toledo GmbH, Laboratory & Weighing Technologies, CH-8606 Greifensee, Switzerland