Quick Setup Guide

InTrac 785/787 Retractable Housings

Contents

cs	Průvodce rychlým nastavením	2
da	Lynvejledning	5
de	Kurzanleitung	8
en	Quick Setup Guide	11
es	Guía de configuración rápida	14
fi	Pika-asetusopas	17
fr	Guide de paramétrage rapide	20
hu	Gyorsbeállítási útmutató	23
it	Guida alla configurazione rapida	26
ja	クイック セットアップ ガイド	29
ko	사용 매뉴얼	32
nl	Beknopte handleiding	35
pl	Podręcznik szybkiej konfiguracji	38
pt	Guia de Configuração Rápida	41
ru	Руководство по быстрой настройке	44
sv	Snabbguide	47
th	คู่มือการตั้งค่าอย่างรวดเร็ว	50
zh	快速设置指南	53

InTrac 785/787

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1 Úvod

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Quick Setup Guide

1 Introduction

The METTER TOLEDO InTrac™ 785/787 retractable housings are intended for use in conjunction with specified METTLER TOLEDO electrodes/sensors, namely pH and redox (ORP) combination electrodes as well as dissolved oxygen (DO), conductivity or turbidity sensors. Please read through this quick set-up guide and the operating instructions carefully before commissioning, in order to ensure trouble-free use. Operating and maintenance should be carried out only by trained personnel and staff who have read and understood the operating instructions.

2 Safety instructions

 **Note:** Before every start-up, the housing must be checked for:

- Damage to the connections, fastenings, etc.
- Leakage
- Perfect functioning
- Authorization for use in conjunction with other plant equipment and resources.

 **Note:** The manufacturer/supplier accepts no responsibility for any damage caused by unauthorized attachments or for the incorporation of spare parts which are not of METTLER TOLEDO provenance. The risk is borne entirely by the operator.

 **Caution:** Before the housing is put into operation, the operator must have already clarified that use of the sensor in conjunction with the other associated equipment and resources is fully authorized.

 **Caution:** A defective housing may neither be installed nor put into operation. Faulty containment or installation out of conformance with regulations and instructions can lead to the escape of medium or to pressure surges (explosion), potentially harmful both to persons and to the environment.

 **Caution:** The housing may only be operated together with the specified electrodes/sensors. If an electrode/sensor is broken it must be replaced immediately.

 **Warning:** Before dismantling the housing or commencing any maintenance work on it, ensure that the equipment in which the housing is installed is in a safe condition (depressurize, no explosion risk, empty, rinse, vent, etc.).

 **Warning:** The electrode or sensor may only be removed from a mounted retractable housing when this is in the "service" (retracted/withdrawn) position and the ball valve is closed, since if the electrode/sensor is missing from the housing there will be path for the process medium to escape to the environment.


 **Warning:** The housing may only be moved to the "measure" position when a sensor is fitted and the ball valve is open. The maximum permissible temperature and pressure specifications must not be exceeded. The values depend upon the design and type of the retractable housing. The relative specifications are given on the label of the individual housings.

3 Product description

The label on the housing contains the following information

METTLER TOLEDO www.mt.com Im Hackacker 15 CH-8902 Urdorf, Switzerland	Manufacturer of the housing and contact details
P/N	Part number
S/N	Serial number
PS	Maximum permissible pressure
TS	Maximum permissible temperature
CE _{xxxx}	CE marking with notified body number
See instruction manual	See instruction manual for detailed description

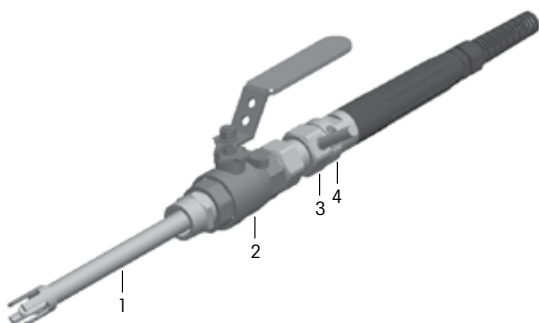4 Installation and commissioning

1. While unpacking, check the housing for mechanical damage. Please immediately report any damage to your local METTLER TOLEDO representative.
2. Ensure the housing is fully retracted and in the "service" position.
3. Isolate the pipe/tank/vessel etc. from the housing (depressurize, empty, rinse, purge and vent, etc.) before starting any installation work.
4. Fit the housing to flange/NPT thread of process connection.
5. Attach the flushing lines (optional flushing chamber when applicable).
6. Fit the electrode/sensor to the housing and connect the electrode/sensor and the transmitter with corresponding cable.
7. Loosen fixing screw **<3>** to release the immersion tube **<1>** by hand or using a no. 38 spanner.
8. Open the ball valve **<2>**.
9. Slide the immersion tube into the measuring position as far as necessary or until it stops.
10. If the housing is not inserted as far as the stop, fix the adjusting ring **<4>** in the desired position.
11. Tighten fixing screw **<3>** by using a no. 38 spanner.

Refer to housing instruction manual for detailed description.

5 Operation: Inspection work

1. Visually check fastenings (ring nut, flange, NPT thread) of the housing at the process connection for firm sealing and leaks.
2. If necessary, check the condition of the electrode/sensor (refer to sensor instruction manual for cleaning and calibration procedures).
3. Calibrate/clean the electrode/sensor if necessary (refer to sensor and transmitter instruction manuals).

6 Maintenance

1. Ensure the housing is fully retracted and in the "service" position.
2. Isolate the pipe/tank/vessel etc. from the housing (depressurize, empty, rinse, purge and vent, etc.).
3. Remove the sensor from the housing.
4. Remove the retractable housing from the pipe/tank/vessel etc. (if changing of housing seal is needed).
5. Change the housing seal.
6. Calibrate/clean the electrode/sensor with reference to sensor and transmitter instruction manuals.
7. Follow the installation and commissioning procedure.

7 Environmental protection

Waste electrical products should not be disposed of with household waste. Please recycle where facilities exist. Check with your Local Authority or retailer for recycling advice.

8 Warranty

On manufacturing defects, 12 months after delivery.

9 Storage condition

We recommend following the shutdown procedure as described in the housing instruction manual. The cleaned and dried housing should be stored at room temperature.

10 EC Declaration of Conformity

The complete declaration is available in the declaration of conformity certificate

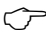
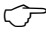
InTrac is a trademark of the METTLER TOLEDO Group.

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All rights reserved