# iLink Multi

## Instruction Manual

cs	Návod k obsluze	2	ko	사용 매뉴얼	42
da	Brugsanvisning	6	nl	Gebruikershandleiding	46
de	Betriebsanleitung	10	pl	Instrukcja obsługi	50
en	Instruction manual	14	pt	Manual de instruções	54
es	Manual de instrucciones	18	ru	Инструкция по эксплуатации	58
fi	Käyttöohje	22	sv	Bruksanvisning	62
fr	Instructions d'utilisation	26	th	คู่มือคำแนะนำ	66
hu	Használati utasítás	30	zh	说明书	70
it	Istruzioni per l'uso	34			
ja	取扱説明書	38			

iLink Multi  
30 370 185 C

**METTLER** **TOLEDO**

## 1 Úvod

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## 1 Introduction

The iLink™ Multi is a universal device for connecting digital ISM® sensors (1-wire; RS485) to a PC/laptop running iSense™ software.

iLink Multi contains precision ambient pressure and relative humidity sensors for simplifying the calibration process of oxygen and dissolved oxygen sensors. Also, a connection terminal to attach an external temperature probe is available to increase calibration convenience.

## 2 Technical data iLink Multi

Measurement parameters	pH, amperometric gas and dissolved oxygen, optical dissolved oxygen (oDO), CO <sub>2</sub>
------------------------	--

Communication to PC	USB (cable included)
---------------------	----------------------

Sensor communication	1-wire (AK9/BNC50 cable not included)* RS485 (iLink Multi cable/Set oDO not included)**
----------------------	--

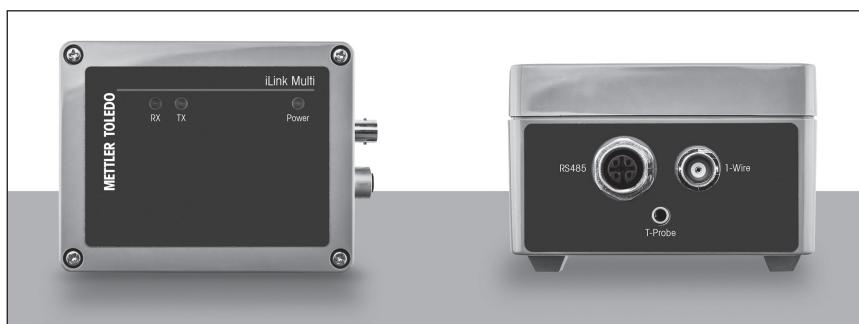
Relative humidity	5 ... 95 % rh (non-condensing)
-------------------	--------------------------------

Environmental temperature	-15 to +50 °C
---------------------------	---------------

\* P/N 59 902 168 to order the AK9 / 1 m/BNC50 cable

\*\* P/N 30 355 582 to order the iLink Multi cable/Set oDO (RS485)

## 3 Operating elements

### 3.1 RS485

① Use the 5-pin to VP8 adapter for VP8 connecting sensors

② Use the 5-pin cable for 5-pin (M12) connecting sensors

### 3.2 1-wire

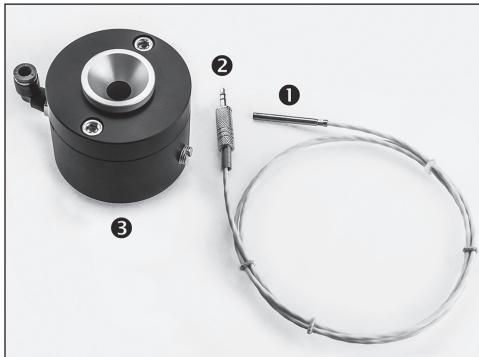
① Connection to iLink Multi (male BNC50)

② Connection to 1-wire sensor

### 3.3 Jack connection to temperature probe

The iLink Multi offers the possibility to connect an external temperature probe. This shortens the reaction time when calibrating optical dissolved oxygen sensors.

When using the METTLER TOLEDO CalBox™ (P/N 52 300 400) an appropriate temperature probe is included.

① Temperature probe

② 2.5 mm Jack connection to iLink Multi

③ CalBox

① CalBox with mounted temperature probe

**Note: Remove protective cap from the CalBox**

The temperature probe must be inserted gently to reach its measurement position.

## 4 Connection to iSense

Use the supplied USB cable to connect the iLink Multi to a PC/laptop.

ISM is a registered trademark of the Mettler-Toledo Group in Switzerland, Brazil, USA, China, European Union, South Korea, Russia and Singapore.

CalBox, iLink and iSense are trademarks of the Mettler-Toledo Group.

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All rights reserved