T100 M12–InTap BT Sensor Interface Instruction Manual

–	Regulatory approval	2
cs	Návod k obsluze	4
da	Brugsanvisning	6
de	Betriebsanleitung	8
en	Instruction manual	10
es	Manual de instrucciones	12
fi	Käyttöohje	14
fr	Instructions d'utilisation	16
hu	Használati utasítás	18
it	Istruzioni per l'uso	20
ja	取扱説明書	22
ko	사용 매뉴얼	24
nl	Gebruikershandleiding	26
pl	Instrukcja obsługi	28
pt	Manual de instruções	30
ru	Инструкция по эксплуатации	32
sv	Bruksanvisning	34
th	คู่มือคำแนะนำ	36
zh	说明书	38

T100 M12–InTap BT Sensor Interface
30 448 013 B

Regulatory approval

FCC Statement

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy, and if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

IC Statement

This device complies with Industry Canada license-exempt RSS standard(s). Operation is subject to the following two conditions: (1) this device may not cause interference, and (2) this device must accept any interference, including interference that may cause undesired operation of the device.

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Regulatory approval

For regulatory approvals please refer to pages 2 to 3.

1 Introduction

The T100 is a Bluetooth interface accessory for the InTap™ portable optical dissolved oxygen analyzer. It can be installed between a transmitter (M400/M800) and a METTLER TOLEDO optical dissolved oxygen sensor using the M12 connector.

Once installed, the InTap analyzer can communicate calibration values through the Bluetooth interface and hence calibrate the attached sensor.

The communication between sensor and transmitter is not interrupted by this installation.

2 Technical data T100

Measurement parameters	RS485: optical DO
Transmission frequency	2.4 GHz, Bluetooth 4.0
Range	10 m
Power supply	Provided by connected transmitter M400/M800 (24 V, min 375 mA)
Operating temperature	–15 to +60 °C
Relative humidity	5 ... 95 % rH (non-condensing)
Protection class	IP67

2.1 Connecting sensors

The T100 connects in between a sensor and an M400/M800 transmitter. The T100 is installed on top of the sensor and its output connects to the transmitter cable.

3 Pairing with InTap

The interface works in conjunction with the Bluetooth-enabled InTap optical dissolved oxygen analyzer.

An in-line sensor that uses a T100 interface can be calibrated wirelessly. Additionally, the InTap stores historical and new calibration date of the installed in-line sensors.

4 LED messages

LED green on	Powered up and connected to sensor
LED blue on	Bluetooth connection established between the InTap and the in-line sensor
LED yellow on	At least one warning registered. See diagnostics in iSense iMonitor.
LED red on	Hardware error or communication interrupted

5 Environmental protection

Waste electrical products should not be disposed of with household waste. Please recycle where facilities exist. Check with your Local Authority or retailer for recycling advice.

InTap and iSense are trademarks of the METTLER TOLEDO Group.

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All rights reserved