METTLER TOLEDO DM141-SC

Silver ring electrode
Silberring-Elektrode
Electrode à l'anneau d'argent
Electrodo con anillo de plata
Электрод с серебряным кольцом
金属银环电极

Combined silver ring electrode
with ceramic frit for argentometric titrations.

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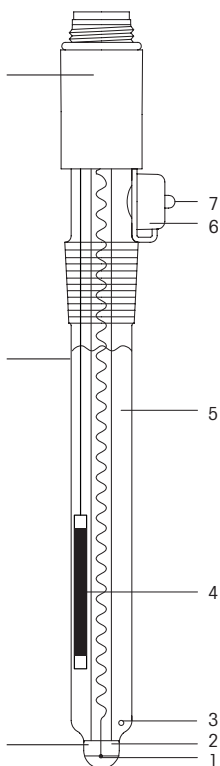
1. Pre-operation treatment

- 1.1 To equalize pressure remove rubber cap (6) or perforate its nipple (7).
- 1.2 Refill the reference electrolyte METTLER TOLEDO No. 51340047 to approx. 1 cm below the aperture (6).
- 1.3 Place the electrode in deion. water for 15 minutes to dissolve the KNO_3 crystals in or on the frit (3).

2. Testing

The electrode can be tested by measurement of its potential with a 0.1 mol/L AgNO_3 test solution and its response time or a titration.

- 2.1 The electrode potential for the 0.1 mol/L AgNO_3 test solution should be within approx. 250 and 350 mV.
- 2.2 Response time of the electrode: After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 2.3 A titer determination of 0.1 mol/L AgNO_3 is best suited for a test titration. The titration has to show a sharp potential jump of at least approx. 130 mV. The equivalence point should be in the range of 50–200 mV.

3. Storage and lifetime

- 3.1 Store the electrode in the reference electrolyte – the frit has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 3.2 Never let the electrode get dry, otherwise the KNO_3 -crystals in or on the frit have to be dissolved.
3. If testing a used electrode the requirements indicated under 2.3 should hold true.

4. Maintenance Instructions

- 4.1 The tolerable temperature range extends from 0 to 80 °C (32–176 °F).
- 4.2 The sample should never enter through the frit into the reference electrode. Therefore the level of the reference electrolyte must always be higher than the sample solution level.
- 4.3 Solutions containing protein quickly block the frit, because protein precipitates on it.
- 4.4 Do not let the cable connector plug corrode.
- 4.5. Ensure that your sample is acidified with diluted nitric or citric acid. Strictly avoid highly concentrated nitric acid, which leads to corrosion of the silver ring!

5. Cleaning procedures

- 5.1 Contamination of the electrolyte: drain out the electrolyte solution and refill with a new one METTLER TOLEDO No. 51340047. Place the electrode for several hours in deion. water.
- 5.2 Blocked frit by protein: place it at least for 1 hour in pepsin solution (5% pepsin in 0.1 mol/L HCl, e.g. Article No. 51340068).
- 5.3 Contamination of the silver surface: depending on the contamination several cleaning procedures can be carried out:
 - a) Place the electrode for a few minutes in an ultrasonic bath with deion. water or ethanol.
 - b) Place the electrode for 60 s in sulfochromic acid or 10% HNO₃ and rinse it thoroughly with deion. water.
 - c) Slightly polish the silver surface with Al₂O₃ powder.

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