METTLER TOLEDO

DGI101-SC/DGI102-Mini

Glass electrode with Sensor Chip
Glaselektrode mit Sensor Chip
Electrode de verre avec Sensor Chip
Electrodo de vidrio con Sensor Chip
Стеклянный электрод
с технологией Sensor Chip
具有智能芯片的玻璃电极

Plug & Play combined pH electrode
with ceramic frit for direct pH measurements
and acid / base titrations in aqueous solu-
tions – for small sample volumes.

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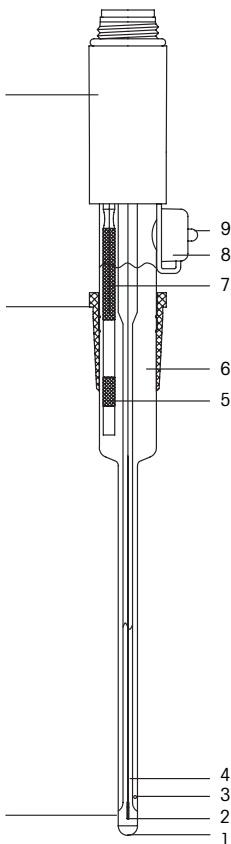
1. Start-up maintenance

- 1.1 The internal buffer solution (4) must fill the entire inside of the glass membrane (1). Remove air bubbles by shaking the electrode gently.
- 1.2 To equalize pressure remove the rubber cap (8) or perforate its nipple (9).
- 1.3 Refill the reference electrolyte (6) – (e.g. ME-51340045) – just up to the aperture (8). Ensure that the silver chloride reference element (7) is always immersed in the reference electrolyte (6).
- 1.4 To activate the glass membrane place the electrode in a tenfold diluted reference electrolyte solution. Use a 0.5 mol/L KCl solution if no reference electrolyte is at hand.

2. Plug & Play performance

The Sensor Chip which stores the sensor parameter is located in the screw cap of the sensor.

The sensor is recognized after connection to the sensor input of the titrator as well as after each restart of the system by reading the sensor parameter on the chip. A popup appears on the titrator terminal when the sensor was successfully recognized. This popup allows you to directly enter the sensor setup. You find further information concerning the functionality of the Plug & Play sensor.

3. Adjustment and calibration

The parameters zero-point and slope are measured by adjustment. A new electrode should have parameter values according to 3.1– 3.2. Instructions for the adjustment method are given in the titrator manual.

- 3.1 The zero point of the electrode (pH value at $E = 0$ mV) is approx. pH 7. The electrode potential, measured in a buffer solution of pH 7, must be between -30 mV and $+30$ mV.
- 3.2 The electrode slope (potential change per pH unit) at the pH range 4–9 must be greater than -55 mV/pH (at 25 °C/ 77 °F).
- 3.3 You should test the response time of the electrode. After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 3.4 If calibrating a used electrode the zero point should be within pH 6 and 8, the slope not less than -52 mV/pH, and the potential change after 60 s should not exceed 3 mV/30 s.

4. Storage and lifetime

- 4.1 Store the electrode in the reference electrolyte – the frit (3) has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 4.2 Never let the glass membrane dry out, otherwise it has to be reactivated and the KCl crystals in/on the frit have to be dissolved.

- 4.3 The lifetime of the electrode is 6 months to 3 years, depending on its use and treatment.

5. Maintenance instructions

- 5.1 The measuring range extends from pH 0 to 14, the tolerable temperature range from 0 to 100 °C (32 to 212 °F).
- 5.2 The sample solution should never enter into the reference electrode through the ceramic frit. Therefore the level of the reference electrolyte must always be higher than the sample solution level.
- 5.3 No reaction should take place between the sample solution and the reference electrolyte: sulfides, bromides, iodides and cyanides react with Ag^+ and some cations (Ag^+ , Hg^{2+} , Au^{3+} , Pb^{2+}) react with Cl^- of the reference electrolyte and cause a blockage of the frit.
- 5.4 Proteinous solutions will quickly block the frit (3) due to protein depositions. To delay this occurrence and therefore to extend the useful measuring time, only use the original reference electrolyte (3 mol/L KCl solution, free of silverchloride).
- 5.5 Just blot the electrode after rinsing, never rub it dry.
- 5.6 Do not let the cable connector plug corrode.
- 5.7 A dried out glass membrane can be reactivated by placing it for several hours in 0.1 mol/L HCl.

6. Cleaning procedures

- 6.1 Contamination of the reference electrolyte: Drain the reference electrolyte through the aperture (8) and proceed according to 1.3 and 1.4.
- 6.2 Frit blocked by AgCl : place it overnight in concentrated ammonia, rinse and condition it for 1 hour in a buffer solution pH 4.
- 6.3 Frit blocked by sulfide precipitate: place it in a solution of thiourea (8% thiourea/1 mol/L HCl, e.g. METTLER TOLEDO No. 51340070) until the dark color has vanished.
- 6.4 Frit blocked by protein: place it at least for 1 hour in pepsin solution (5% pepsin/0.1 mol/L HCl, e.g. METTLER TOLEDO No. 51340068).
- 6.5 Other contaminations: place it for a few minutes in an ultrasonic bath with deion. water, ethanol or a mixed acid solution (e.g. aqua regia).

7. Regeneration

If the electrode does not meet the qualifications according to 3.4 anymore, try a corrosion of the glass membrane: place it 1–2 minutes in diluted HF-solution (2% HF/5% HCl, e.g. ME-51340073), rinse thoroughly with deion. water and treat it as in 6.1.

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