METTLER TOLEDO

DM1102-SC

Silver ring electrode with Sensor Chip
Silberringelektrode mit Sensor Chip
Electrode combinée à anneau d'argent
avec Sensor Chip
Electrodo con anillo de plata combinado
con Sensor Chip
Электрод с серебряным кольцом
с технологией Sensor Chip
具有智能芯片的金属银环电极

Plug & Play combined silver ring electrode
 with ceramic frit for argentometric titrations –
 for small sample volumes.

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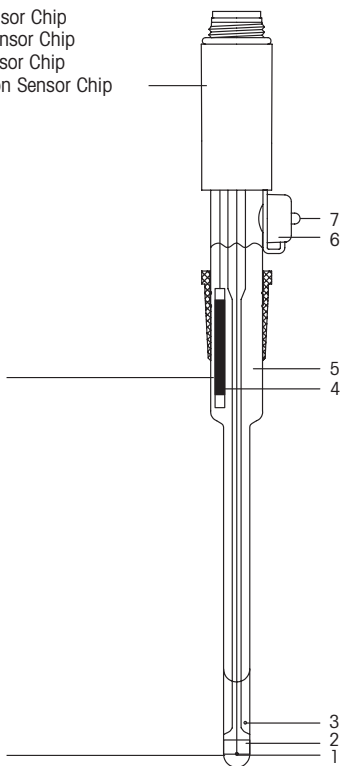
1. Start-up maintenance

- 1.1 To equalize pressure remove rubber cap (6) or perforate its nipple (7).
- 1.2 Refill the reference electrolyte METTLER TOLEDO No. 51340047 to approx. 1 cm below the aperture (6).
- 1.3 Place the electrode in deion. water for 15 minutes to dissolve the KNO_3 crystals in or on the frit (3).

2. Plug & Play performance

The Sensor Chip which stores the sensor parameter is located in the screw cap of the sensor.

The sensor is recognized after connection to the sensor input of the titrator as well as after each restart of the system by reading the sensor parameter on the chip. A popup appears on the titrator terminal when the sensor was successfully recognized. This popup allows you to directly enter the sensor setup. You find further information concerning the functionality of the Plug & Play sensor.

3. Testing

The electrode can be tested by measurement of its potential with a 0.1 mol/L AgNO_3 and its response time or a titration.

- 3.1 The electrode potential for the 0.1 mol/L AgNO_3 test solution should be within approx. 250 and 350 mV.
- 3.2 Response time of the electrode: After 30 s the indicated electrode potential should not change more than 2 mV within the following 30 s.
- 3.3 A titer determination of 0.1 mol/L AgNO_3 is best suited for a test titration. The titration has to show a sharp potential jump of at least approx. 130 mV. The equivalence point should be in the range of 50–200mV.

4. Storage and lifetime

- 4.1 Store the electrode in the reference electrolyte – the frit has to be immersed – and close the aperture (perforated rubber cap is sufficient).
- 4.2 Never let the electrode get dry, otherwise the KNO_3 crystals in or on the frit have to be dissolved.
- 4.3 If testing a used electrode the requirements indicated under 3.3 should hold true.

5. Maintenance Instructions

- 5.1 The tolerable temperature range extends from 0 to 80 °C (32 – 176 °F).
- 5.2 The sample should never enter through the frit into the reference electrode. Therefore the level of the reference electrolyte must always be higher than the sample solution level.
- 5.3 Solutions containing protein quickly block the frit, because protein precipitates on it.
- 5.4 Do not let the cable connector plug corrode.
- 5.5 Ensure that your sample is acidified with diluted nitric or citric acid. Strictly avoid highly concentrated nitric acid, which leads to corrosion of the silver ring!

6. Cleaning procedures

- 6.1 Contamination of the electrolyte: drain out the electrolyte solution and refill with a new one METTLER TOLEDO No. 51340047. Place the electrode for several hours in deion. water.
- 6.2 Blocked frit by protein: place it at least for 1 hour in pepsin solution (5% pepsin in 0.1 mol/L HCl, e.g. Article No. 51340068).
- 6.3 Contamination of the silver surface: depending on the contamination several cleaning procedures can be carried out:
 - a) Place the electrode for a few minutes in an ultrasonic bath with deion. water or ethanol.
 - b) Place the electrode for 60 s in sulfochromic acid or 10% HNO₃ and rinse it thoroughly with deion. water.
 - c) Slightly polish the silver surface with Al₂O₃ powder.

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