## Changing the interior body Ersetzen des Innenkörpers Remplacement du corps interne

### InPro® 6800/6900/6950 Series O<sub>2</sub> Sensors 12/25 mm

## Changing the interior body

**Warning!** The O<sub>2</sub> electrolyte has an alkaline pH value of 13. Contact of electrolyte with mucous membrane or eyes is to be avoided. **Therefore protective gloves and safety glasses have to be worn for the following dismantling works.**

If such contact occurs, the affected area should be well rinsed with water. In the case of accident, or should ever any adverse signs appear, get immediate medical attention.

When changing the interior body, please observe the following instructions (see also the following illustration):


**Attention!** Make sure that this maintenance step is carried out in **clean place**.

1. Unscrew the cap sleeve from the sensor shaft and carefully pull it off the sensor.
2. Pull off the membrane body from the interior body. If it is tight-fitted, eject by pushing it with the flat finger tip. Before electrolyte is refilled, the membrane body must be removed from the cap sleeve.
3. Dry the interior body with a paper tissue.
4. Unscrew the retainer nut of the interior body with an adjustable wrench or with a  $\frac{3}{8}$ " or 9 mm (InPro 6850i) wrench.
5. Remove the interior body by pulling it out of the sensor shaft. If necessary use a plier.

**Warning!** Do not twist the interior body. Otherwise the connection pins can be damaged.

6. Insert the new interior body in the sensor shaft. Turn the interior body in the shaft until the slit of the interior body is aligned with the pin placed in the shaft.
7. Press the body in the shaft and screw the new retainer nut in place.
8. Examine the O-rings visually for mechanical defects, and replace if necessary.
9. Half-fill the new membrane body with O<sub>2</sub> electrolyte and make sure that all bubbles are removed. Air bubbles can be removed by carefully tapping on the membrane body.
10. Slip the membrane body over the interior body while holding the sensor in a vertical position. The excess electrolyte will be displaced and have to be removed with a paper tissue.  
**Important!** No electrolyte, sample media or contamination may be present between the membrane body and the cap sleeve. Please check carefully.
11. Carefully slip the cap sleeve over the fitted membrane body, holding the sensor in a vertical position and screw it tight. The cap sleeve must be clean and dry.
12. After each exchange of electrolyte or membrane body, the sensor has to be repolarized and recalibrated.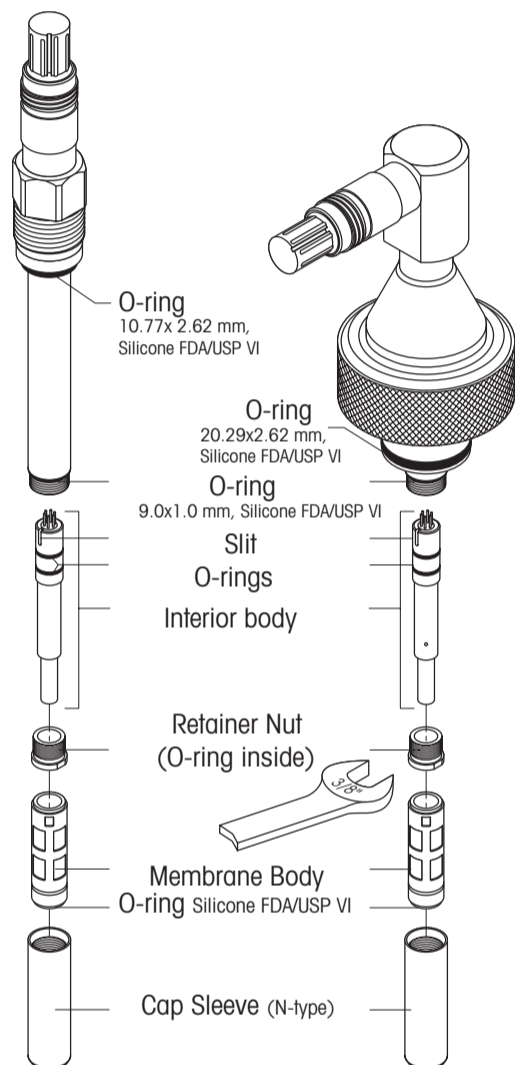

**BR** Mettler-Toledo Ind. e Com. Ltda., Alameda Araguaia, 451, Alphaville, BR-06455-000 Barueri/SP, Brazil, Phone +55 11 4166 7444, Fax +55 11 4166 7401

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