Adjustment for Verification

JL-GE, PL-E

▶ www.mt.com/Adjust-JLGE-PLE-INI

English

Introduction

State or local laws and regulations may require this instrument to be inspected and certified by an authorized inspector or agent before you may use it in commerce with customers.

METTLER TOLEDO takes reasonable steps to keep its products in conformity with various nationally recognized standards and regulations. However, our prices do not include the cost of any related inspections, permits or inspection fees that may be required by any federal, provincial, state or local laws, ordinances, regulations, codes or standards.


Initial adjustment for verification

After putting the balance into operation, an adjustment dialog appears on the display. After executing this adjustment, the dialog disappears.


- The adjustment weight is available.
 - The weighing pan is empty.
- 1 Connect the balance to the power supply using the appropriate AC/DC adapter and country-specific power cable.
 - 2 Level the balance.
 - 3 Press to switch on the balance.
 - 4 Let the balance warm up for approximately 30 minutes.
 - ➔ The display shows the required adjustment weight.

METTLER **TOLEDO**

- 5 If the displayed value does not match your adjustment weight, press Col until the displayed weight matches your adjustment weight.
- 6 Press \leftarrow to confirm (to abort, press $\rightarrow 0 \leftarrow$).
 - The display flashes.
- 7 Place the adjustment weight in the center of the weighing pan.
 - The adjustment process starts.
 - The display flashes **0.00 g**.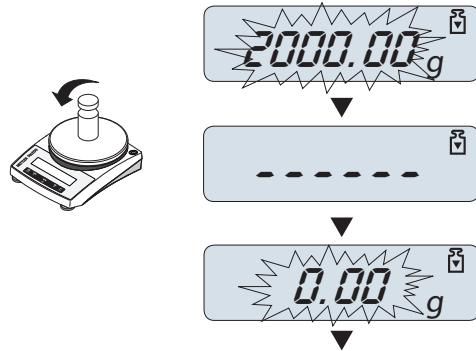

- 8 Remove the adjustment weight.
- When the adjustment is finished, the balance briefly shows **ADJ.DONE**.
- The balance is ready for operation.

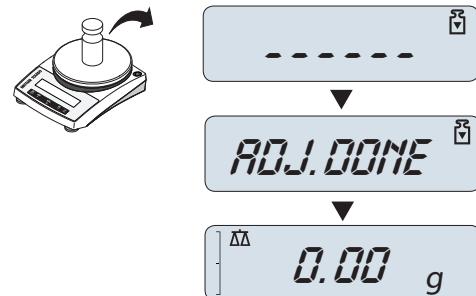Re-adjustment for verification

- 1 Perform steps 1 to 4 of the initial adjustment.
- 2 Press and hold Off to switch off the balance.
- 3 Press the service button (1) on the back of the balance (e.g. using a paperclip) until the display shows **CAL**.
 - The display shows the required adjustment weight.
- 4 Continue from step 5 of the initial adjustment.

Approved balances

After adjustment, the display shows the last digit in brackets (approved-balance mode).

Contact your local weights and measures office to have your balance certified as an approved balance and to have a security seal affixed.

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© 05/2022 METTLER TOLEDO. All rights reserved.
Mettler-Toledo GmbH